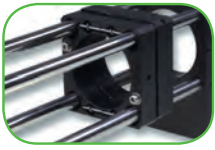
LINOS CATALOG

Edition λ

An Excelitas Technologies Company

Mechanics

High-precision opto-mechanics for reliable integration of optics

Mikrobench
11

Nanobench
79

Tube System C
101

Positioning Systems
127

Profile and Rail Systems
167

Mounts, Posts and Holders
189

Mirror Mounts
223

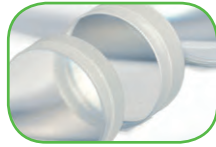
Q-Sets
237

Optics

Optics with superior performance

Singlets
251


Achromats and Lens Systems
283


Plano Optics
295

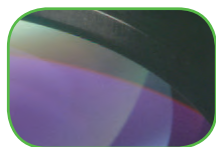Mirrors
331

Thin Film Coatings
359

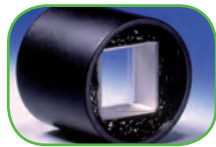Polarization Optics
389

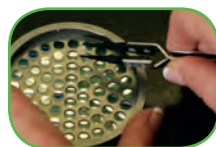
Laser Optics
405

Machine Vision Solutions
423

Micro Imaging Solutions
449

Capabilities
473

Optics Software

Optic Design Software and lens configuration

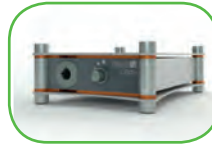
WinLens™ and MachVis
485

Instruments

High-quality laser and light sources for demanding applications

Laser
559

Light Sources
579

Electro Optics

High-performance Faraday isolators, Pockels cells and Modulators

Faraday Isolators
505

Laser Modulators
and Pockels Cells
529

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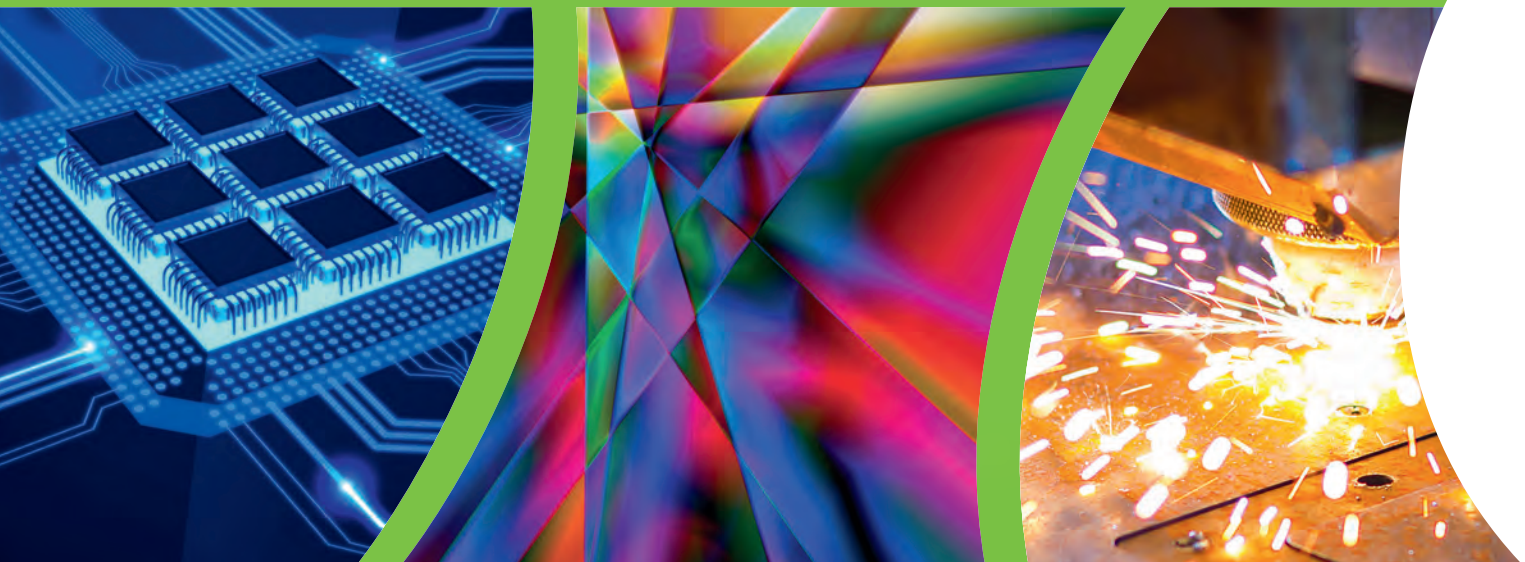
LINOS - Premium brand quality criteria

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Isabel Haackert, Sales Director Catalog & Technical Sales

Dear Readers,

With the new edition λ of the LINOS Catalog, we offer you our extensive range of high-quality standard products. Specifications and useful additional information will help you find solutions for your scientific research requirements, image processing, laser material processing, laser sources and LED illumination and optical measurement technology.

In addition, as a manufacturer of high-end optics we also offer you customized solutions. With our extensive optical fabrication capabilities and design expertise (see page 473), we offer a wide range of optics, from special optics to micro-optics, from aspheres to free form surfaces.

*Use our expertise to meet your requirements! We look forward to receiving your inquiries and being of assistance.
Best regards,*

Isabel Haackert, Sales Director Catalog & Technical Sales

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Excelitas Technologies Corp. is a global technology leader focused on providing innovative, high-performance photonic solutions. Whether you seek off-the-shelf photonic products, or complex custom components, modules and systems, our photonic technologies represent the highest caliber of quality, reliability and performance.

With the acquisition of Qioptiq in 2013, we possess a distinctive brand heritage encompassing pedigrees such as LINOS, Rodenstock, Spindler & Hoyer, Gsanger, Optem, X-Cite, Point Source and OmniCure, amongst others. Our collective knowledge and expertise delivers sophisticated illumination, optical, electro-optical and detection solutions to global customers across the biomedical, scientific research, safety and security, consumer products, semiconductor, industrial manufacturing, and defense and aerospace sectors. Our 5,500 employees operate from 21 state-of-the-art photonics facilities based in 15 different countries across North America, Europe and Asia, and stand committed to fully enabling our customers' innovation and success.

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Visit our online Q-Shop and stay up-to-date! At qioptiq-shop.com you will find up-to-date information about new products, prices and availability. All orders for stock items from Germany that are received working days by 1:00 pm will be sent to you on the same day! You can also request quotes online.

In the Q-Shop you will also find our trade fair and event dates as well as extensive information and numerous publications for download!

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The dialog with our customers is the key to success. Qioptiq employees across three continents are very happy to support you with their expertise and knowledge.

Your Qioptiq contacts

Qioptiq Europe

Phone: +49 (0) 551 6935-0
sales.qioptiq@excelitas.com

Qioptiq North America

Phone: +1 800 775 6786
sales.qioptiq@excelitas.com

Qioptiq Asia Pacific

Phone: +65 6499 7777
sales.qioptiq@excelitas.com

Japan

Opto Science Inc.
Phone: +81 3 3356 1064
info@optoscience.com

Netherlands

VADENO Nederland
Phone: +31 610 414 067
info@vadeno.nl

Mechanics

High-precision opto-mechanics for reliable
integration of optics

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Our quality criteria:

- German manufacturing expertise using only high-quality, extremely durable materials
- Surface finishing with precision anodizing
- Narrow production tolerances
- On-going system expansion with new optical and mechanical components
- The Qioptiq guarantee: Complete compatibility of all Microbench series components - from 1963 to today

LINOS Microbench

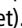
It is our pleasure presenting You the original Microbench, invented by Qioptiq!

This compact 4-rod construction system is in serial production in Göttingen, Germany since 1963. The LINOS Microbench was developed by LINOS predecessor Spindler & Hoyer as an innovative replacement of the triangular rail system. Small in its dimensions, precisely centered, simple to integrate, easy to expand, and very rugged: the LINOS Microbench is ideal for all spatial structures in any compact form.

Thanks to the broad range of optical and mechanical components available, the LINOS Microbench is an integral part of many optomechanical instruments. Take advantage of more than 50 years of experience to ensure the highest precision and stability for your requirements.

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Special Features:

For a rapid setup of your optics, Qioptiq offers a large array of mounted optics developed specially for the Microbench; look for the  symbol in the catalog.

Ideal Areas:

Optics with Ø of 18 to 31.5 mm; illumination and imaging assemblies; optical inspection, measurement and testing equipment; interferometric applications, and more.

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Content

<i>Technical Features</i>	16
Sets	
<i>Microbench Set Mechanics</i>	19
<i>Microbench Set Optics</i>	21
<i>Microbench Basic Set</i>	22
<i>Microbench Set O-UV</i>	24
<i>Laserset</i>	25
<i>Storage Cases</i>	25
<i>Microbench Mechanics Test Kit</i>	26
Mounting Plates	
<i>Mounting Plates 25/30/35/1 "</i>	27
<i>Mounting Plate 25 with Mounting Holes</i>	28
<i>Mounting Plate 25/30/35, Latchable</i>	28
<i>Mounting Plate 25/30, Latchable, 10°</i>	29
<i>Mounting Plate 25 - T10</i>	30
<i>Mounting Plate 25E/30E</i>	30
<i>Mounting Plate M35x0.5, also Latchable</i>	31
<i>Magnetic Cage Assembly Swapper Plates</i>	31
Holder	
<i>Holder 25/30/35/M35x0.5</i>	32
<i>Adjustable Holder 25</i>	32
<i>Rod Holder G/F/90°/V</i>	33
<i>Mounting Bases</i>	34
<i>Rotating Holder, 5 Positions</i>	34
Rods and Cubes	
<i>Rods</i>	35
<i>Cube</i>	35
<i>Beam-Deflector Cube</i>	36
Corner Connectors and Angular Adjustments	
<i>Corner Connector 40/80/40 with Threads</i>	36
<i>Corner Connector T</i>	37
<i>Connector 80 V</i>	37
<i>Swivel Connector</i>	38
<i>Swivel Arm with Scale</i>	38
<i>Mounting Bracket 40</i>	39
Positioners	
<i>XYZ-Fine Adjustment Unit</i>	39
<i>Slip Plate Positioner</i>	40
<i>Centering Mounting Plate/Holder 25</i>	40
<i>X-Y Precision Translation Stage</i>	41
<i>fle.X-Plate</i>	42
<i>Z-Adjustment M</i>	43
<i>Z-Fine Adjustment M</i>	44
<i>Linear Stages for Microbench</i>	45
<i>Z-Translation Stage with Micrometer</i>	46
<i>Z-Translation Stage with Fine-Adjustment Screw</i>	46
<i>Rotary Support</i>	47
<i>Rotary Support 20°</i>	47
<i>Motorized Rotation Mount</i>	48
<i>Counter Support</i>	49
<i>Holder for Micrometer</i>	49
<i>Compression Spring, 60x8x0.7</i>	50

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Content

Fine Adjustment Units and Object Slides

Adjustable Prism Insert 30	50
Adjustable Mirror Insert 30	51
Tilting Platform	51
Prism Table 25, Adjustable	52
Mounting Tool for Beamsplitter Cubes	52
Rotation Adjustment for Cube Inserts	53
Positioning Ring 30	53
Beam Steering Mirror Assembly	54
Optic Support 25	54
Clamp Holder for Prism	55
Prism Support 25	55
Prism Table 25	56
Adjustable Optic Holder 31.5	56

Adapter

Overview Table Adapter	57
Microbench - Nanobench Adapter	58
Adapter 16/25	58
Adapter Collar	59
Cap 30/25	59
Adapter 30/25	59
Adapter 35/30	60
Adapter M35x0.5/30	60
Adapter 35/30-3.5°	60
Adapter 30/C-Mount	61
Adapter 25/C-Mount	61
C-Mount-Adapter-M17	61
Adapter C-Mount/Microbench	62
Adapter 1" to W0.8" or Ø 25 mm	62
Adapter Bench Systems 30 & 1.035"-40 UN	63
Adapter 1.035"-40UN / 30	63
Set of Adapter	64
Adapter 16	64
Adapter 19	64
Adapter 23.2	65
Adapter W0.8"x1/36	65
Adapter W0.8" to 1"	65
Adapter 23.2 to Ø 25	66
Eyepiece Adapter 23.2/30	66
Adapter 25/15	66
FC/FSMA Fiber Adapter	67
FC Fiber Adapter 3.5°	67
Instrument Flange	68
Adapter Plate	68
Set of Retaining Rings	69
Set of Collars	69

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Content

Experiment Set	
Two Beam Interferometer	70
Accessories	
Collimator Holder	71
Cuvette Holder	71
Mounting Plate FLS	72
Closing Caps	72
Focusing Collimators	73
Tube	74
Lens Turret 4x with C-Mount Adapter	74
Microbench Alignment Plate	75
Tools and Screws	
Circlip Kit	75
Special Screwdriver	75
Screwdriver	76
Set of Threaded Pins M2.3x3	76
Set of Threaded Pins M2.3x6	77
Set of Thumbscrews M2.3x9	77
Set of Thumbscrews M2.3x4	77
Set of Screws M2.3x5	77
Set of Hexnuts M2.3	78
Polyamide Plugs	78

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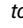
Technical Features

Microbench is a compact component system for numerous areas of application – from lighting and imaging structures to optical experimental setups, measurement / inspection equipment and interferometric applications, to name just a few. The basic principle lies in the combination of standardized mounting plates for integrating optics of all types, connected by stainless steel rods. Both linear and spatial setups can be implemented quickly and easily. All components of the Tube Mounting System C with outer diameter of 30 respectively 35 mm can easily be integrated in the appropriate Microbench mounting plates. Via adapters the LINOS Microbench is also compatible with our Nanobench system.


Qioptiq offers mounted optics developed specially for the Microbench with diameters from 18 mm to 31.5 mm. The special  symbol shows you at a glance which optics are available for your Microbench system.

- Construction system for
 - illumination or imaging systems
 - optical testing, metrology, or inspection systems
 - interferometric systems
 - Compact, three-dimensional setups
- Rigid 4-rod structures
- Optional 2-rod structures for more rapid setup
- Easy to handle
- High mechanical stability

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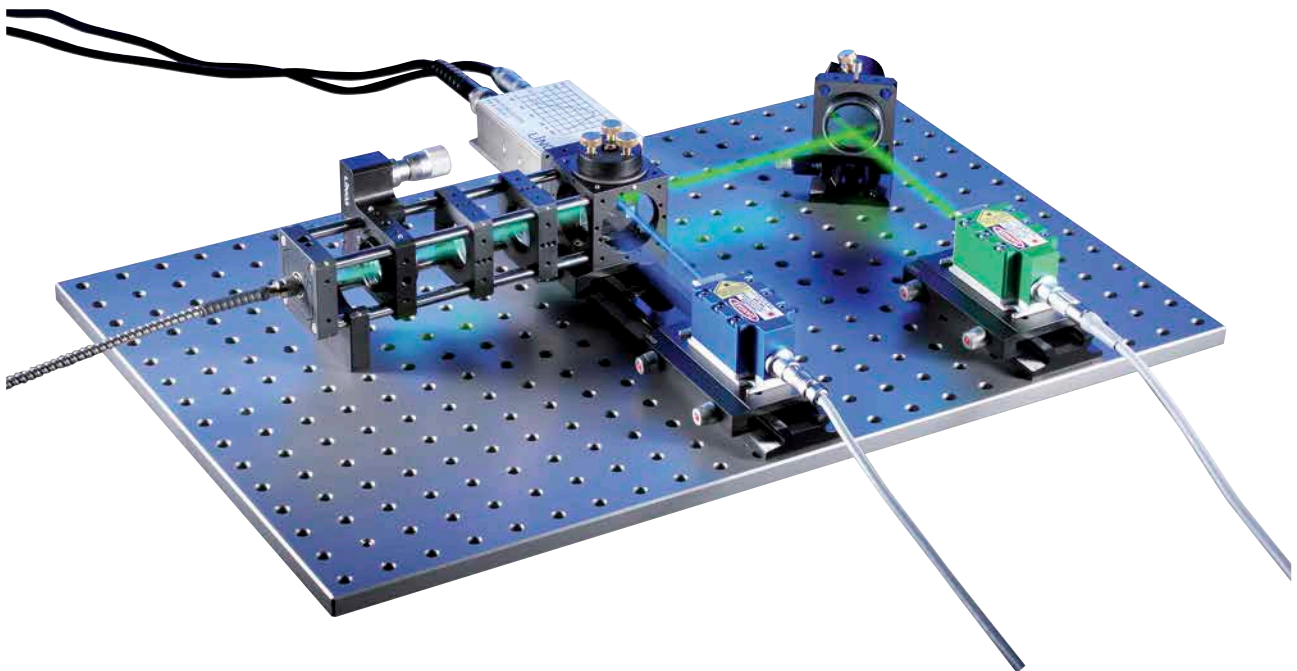Mounting plates

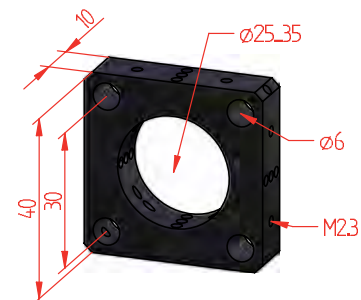
- Lead-free aluminum alloy
- Matt black anodized surfaces
- Inner diameter: 25, 30 or 35 mm with H7 tolerances
- Four bore holes for mounting the connecting rods. The tolerances for the rod bore holes on the mounting plates, together with the tolerances for the rods, guarantee a system with virtually zero play.

Rods


- Hardened stainless steel with polished surfaces
- Lapped to fit
- Outer diameter: 6 mm
- Lengths: 20 mm to 450 mm
- Compatible adjustment rings as holders or stops; e.g., for Z-adjustment of individual components

Cubes

- Multifunctional elements for 3D structures
- Ideal for beam deflectors
- Inner diameter of bore holes: 25 or 30 mm
- Two of the cube sides have bore holes for rods, four sides have threads for mounting plates with tapped holes

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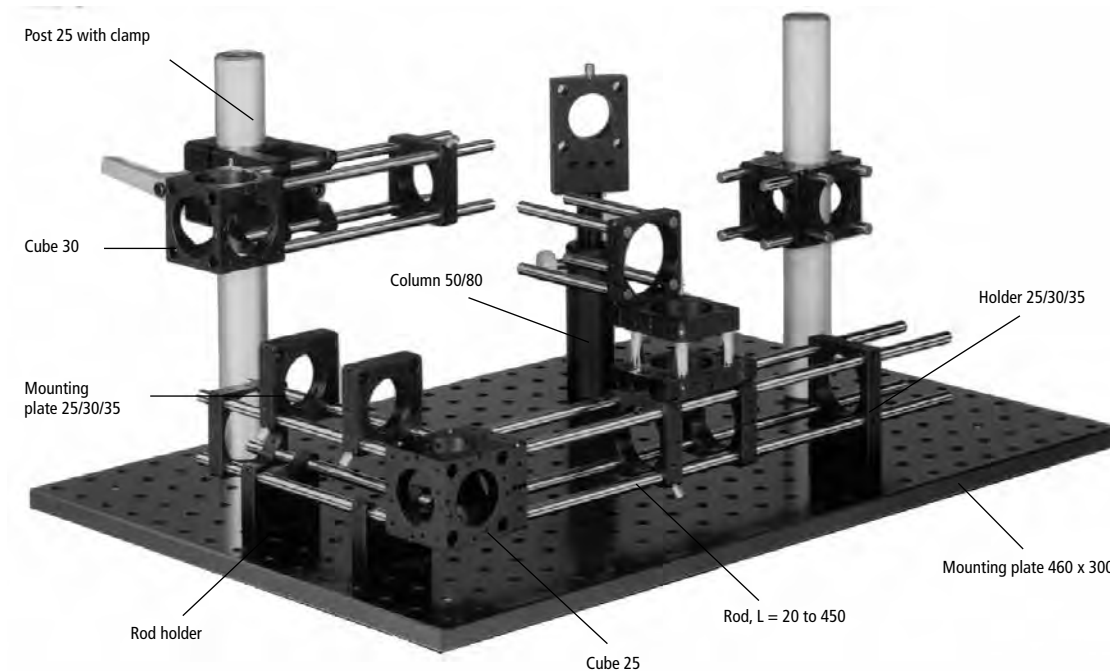- Includes all most important parts for an easy start with Microbench system
- Complete with rails and carriers of the Rail System FLS 40
- Contains tools and installation material

Microbench Set Mechanics

Part No.	Quantity	Description
G06 1355 000	1	Rail FLS 40-250
G06 1374 000	1	Carrier FLR 40-40
G06 1372 000	1	Carrier FLR 40-20
G06 1371 000	2	Carrier FLR 40-10
G06 1309 000	1	Mounting Plate 220x40 mm
G06 5061 000	2	Holder 25
G06 1225 000	2	Holder 30
G06 5084 000	2	Holder 35
G06 1010 000	2	Mounting Plate 25
G06 1042 000	3	Mounting Plate 30
G06 1047 000	2	Mounting Plate 35
G06 1020 000	1	Mounting Plate 25, latchable
G06 1021 000	1	Mounting Plate 30, latchable
G06 1041 000	2	Mounting Plate with mounting holes
G06 1025 000	2	Centering Mounting Plate 25
G06 1080 000	1	Cube 25
G06 1081 000	1	Cube 30
G06 1228 000	1	Rod Holder
G06 1229 000	1	Rod Holder G
G06 1227 000	2	Rod Holder F
G06 1110 000	4	Corner Connector 40 mm
G06 1161 000	1	Counter Support
G06 1160 000	1	Holder for micrometer G 10
G06 1162 000	1	Micrometer G 10
G06 1159 000	2	Compression Spring
G06 5058 000	1	Rotary Support
G06 3730 000	1	Adjustable Mirror Insert 30
G06 3731 000	1	Adjustable Prism Insert 30
G06 5002 000	1	Optic Support 25
G06 1661 000	1	Adapter 1" to W 0.8 and ø 25 mm
G06 1601 000	1	Cap 30/25
G06 1621 000	1	Adapter W0.8x1/36"
G06 1636 000	1	Adapter 23.2
G06 1602 000	1	Adapter 30/25
G06 1658 000	1	Adapter 30/C-Mount
G06 1637 000	1	Adapter 35/30
G06 1650 000	1	Iris Diaphragm C 18 with tube 25
G06 1220 000	1	Set of 10 collars
G06 5092 000	1	Positioning Ring 30
G06 1207 000	4	Rod, 20 mm
G06 1208 000	4	Rod, 30 mm
G06 1209 000	4	Rod, 75 mm
G06 1210 000	4	Rod, 150 mm
G06 1205 000	4	Rod, 200 mm

Microbench Set Mechanics

Item Title	Part No.
Microbench Set Mechanics	G060004000

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- Fast beginning of experiments ensured by wide range of mounted optics

Microbench Set Optics

Part No.	Quantity	Description
G06 3020 000	1	Biconvex lens 20/21.4 mounted
G06 3024 000	1	Biconvex lens 50/21.4, mounted
G06 3026 000	1	Biconvex lens 80/21.4, mounted
G06 3028 000	1	Biconvex lens 150/21.4, mounted
G06 3852 000	1	Biconvex lens 50/24, mounted
G06 3854 000	1	Biconvex lens 100/24, mounted
G06 3014 000	1	Biconvex lens 70/30, in mount 31.5
G06 3042 000	1	Plano-convex lens 30/17, mounted
G06 3043 000	1	Plano-convex lens 40/21.4, mounted
G06 3044 000	1	Plano-convex lens 50/21.4, mounted
G06 3045 000	1	Plano-convex lens 60/21.4, mounted
G06 3046 000	1	Plano-convex lens 80/21.4, mounted
G06 3047 000	1	Plano-convex lens 100/21.4, mounted
G06 3048 000	1	Plano-convex lens 150/21.4, mounted
G06 3049 000	1	Plano-convex lens 200/21.4, mounted
G06 3821 000	1	Plano-convex lens 50/24, mounted
G06 3822 000	1	Plano-convex lens 80/24, mounted
G06 3823 000	1	Plano-convex lens 100/24, mounted
G06 3826 000	1	Plano-convex lens 300/24, mounted
G06 3052 000	1	Plano-convex lens 100/30, in mount 31.5
G06 3057 000	1	Biconcave lens -20/21.4, mounted
G06 3060 000	1	Biconcave lens -40/21.4, mounted
G06 3062 000	1	Biconcave lens -100/21.4, mounted
G06 3892 000	1	Biconcave lens -50/24, mounted
G06 3070 000	1	Plano-concave lens -40/21.4, mounted
G06 3871 000	1	Plano-concave lens -50/24, mounted
G06 3072 000	1	Plano-concave lens -100/21.4, mounted
G06 3076 000	1	Concave-convex lens 50/21.4, mounted
G06 3127 000	1	Achromat 40/17, mounted
G06 3141 000	1	Achromat 50/21.4, mounted
G06 3212 000	1	Achromat 60/24, in mount 25.4
G06 3143 000	1	Achromat 80/21.4, mounted
G06 3144 000	1	Achromat 100/21.4, mounted
G06 3215 000	1	Achromat 120/24, in mount 25.4
G06 3204 000	1	Achromat 160/30, in mount 31.5
G06 3148 000	1	Achromat 200/21.4, mounted
G06 3191 000	1	Achromat -50/17, mounted
G06 3010 000	1	Condenser combination 16/21.4
G06 3099 000	1	Asph. condenser lens 18/21.4, mounted
G06 3098 000	1	Asph. condenser lens 27/30 in mount
G06 3450 000	1	Colored glass filter BG 38, mounted
G06 3451 000	1	Colored glass filter BG 18, mounted
G06 3082 000	1	Colored glass filter VG 9, mounted
G06 3081 000	1	Colored glass filter RG 630, mounted
G06 3080 000	1	Heat absorbing filter KG 1, mounted
G06 3410 000	2	Polarization filter
G06 3411 000	1	Quarter-waveplate
G06 3619 000	1	Front surface concave mirror 14/30 mounted
G06 3628 000	1	Front surface plane mirror ø 30 mounted
G34 0060 000	1	Front surface plane mirror 20 x 30 mm
G34 4143 000	1	Beamsplitter plate 20 x 30 mm
G39 0137 000	1	Plane plate 20 x 30 x 2.5 mm
G33 9913 000	1	90°-Prism 20 x 20 mm
G33 5520 000	1	Beam splitter cube 20 x 20 x 20 mm
G03 8771 000	1	Microscope objective 10x/0.25

Microbench Set Optics

Item Title	Part No.
Microbench Set Optics	G062004000

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Microbench Basic Set

The case contains all components of the following sets:

Microbench Set Mechanics G060004000
Microbench Set Optics G062004000

- Types of systems that may be assembled:
 - Microscopes
 - Off-axis viewers
 - Telescopic systems
 - Measuring telescopes
 - Collimators
 - Polarizers
 - Projection systems
 - Graticule projectors
 - Beamsteering systems
 - Beamsplitter systems
 - Interferometers
 - Illumination systems
- Included: Software Package WinLens™ (see chapter Optics Software for further information)

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Please find further information about application examples, guidelines, frequently asked questions and more in our Brochure "Microbench, Bench and Rail systems - Introduction, solutions and applications".

You will find the brochure for download at each Microbench product under "Docs + Drawings".

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
Microbench Set O-UV

Item Title	Part No.
Microbench Set O-UV	G06202000

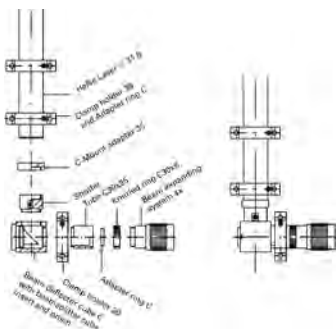Mounting example:
Horizontal angle assembly with beam expander system

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Storage Cases

All storage cases complete with inserts, but without components.

- Suitable for careful and organized storage of Microbench components and components with similar dimensions
- Layout of the compartments as shown for the sets with components

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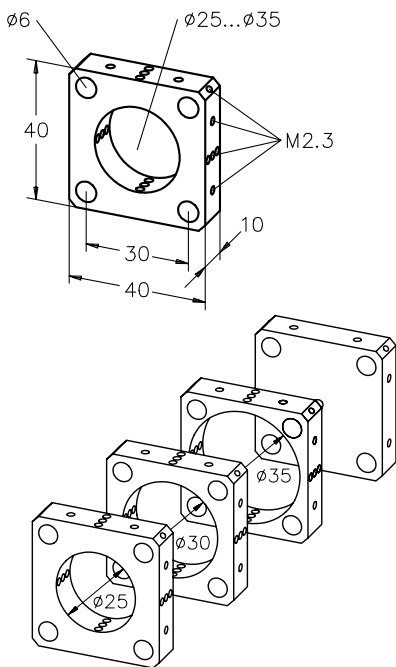
Mounting Plates 25/30/35/1"

Mounting plates are assembled on rods and enable concatenation of multiple optical construction elements. Elements can be positioned as desired along the rod. The threaded pins that secure an element in position exert radial pressure on the rod so the plate position is not altered when the pins are tightened.

All plate elements are made of corrosion-resistant, black anodized, RoHS-compliant aluminum alloy.

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- Mounting plates and holders are delivered without screws. The compatible set of screws are M2.3x3 (qty. 150), part G06 1011 000.
- See section "Fasteners and Tools" for more information on screws

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
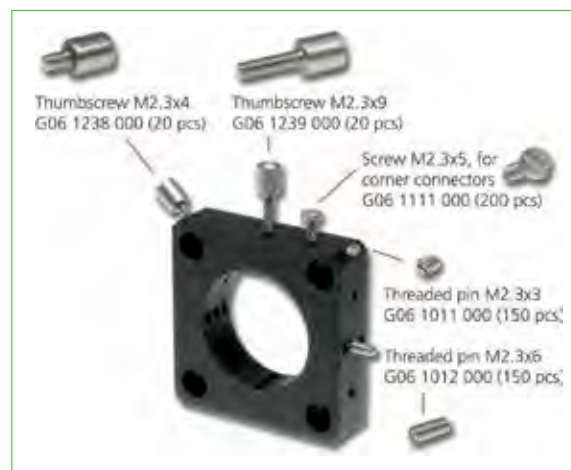
For mounting of optics and components we provide a whole range of screws and threaded pins, please look at section "Fasteners and Tools" and the end of this chapter and add these to your order.

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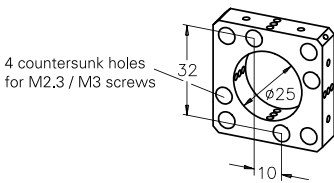Mounting Plate 25 with Mounting Holes



- For attaching to cube 25 or cube 30, for mounting additional rods or attaching to surfaces
- Including screws

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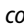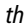For compatible and appropriate optics (lenses, filters, etc), see chapters "Optics". Optics compatible to our bench systems are marked as follows:  for Microbench and  for the Nanobench.

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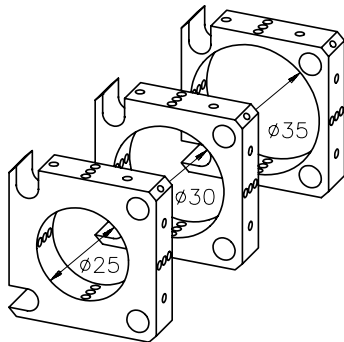
Mounting Plate 25/30/35, Latchable

In contrast to the basic Mounting plates, the latchable models can be temporarily swiveled out from the optical construction, or easily added to an existing construction. The plates are fixed in position with threaded pins or can be held in place by collars.

- Mounting plates are delivered without screws, feel free to order: e.g. Set of screws M2.3x3 (qty. 150) G06 1011 000.
- See section "Fasteners and Tools" for more accessories

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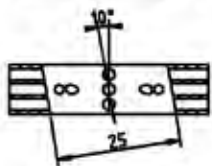
This latchable mounting plate allows the integration of optical components at an 10° angle to prevent back-reflections in the optical path and to prevent etalon effects.

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
- *Mounting plates and holders are delivered without screws. The compatible set of screws are M2.3x3 (qty. 150), part G06 1011 000.*
- *See section "Fasteners and Tools" for more information on screws*

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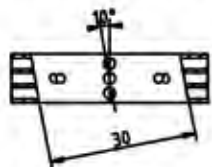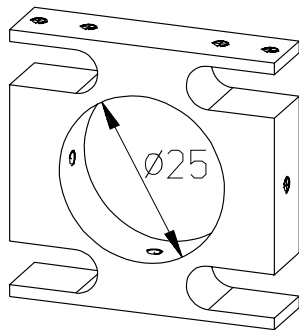
Mounting Plate 25 - T10

- Allows ± 5 mm parallel adjustments when used in combination with T10 / TS10 translation stages
- Please look at the "Positioning Systems" chapter regarding Translation stages T10 / TS10.

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Produktbezeichnung	Part No.
Item Title	Part No.
Mounting Plate 25 - T10	G061070000

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
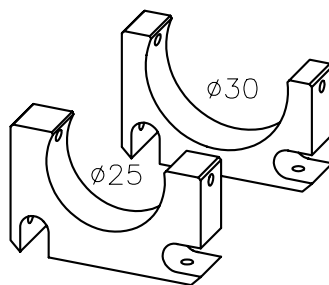
Mounting Plate 25E/30E

- For insertion into existing set-ups

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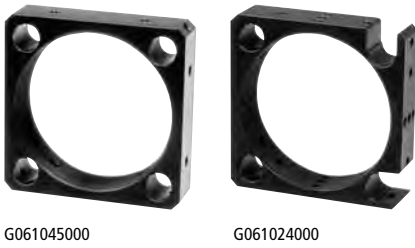G061045000

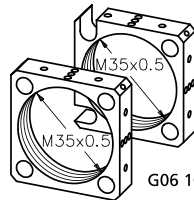
G061024000

Mounting plates M35x0.5, here complete with optics and retaining rings (both not included)

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

G06 1045 000

Mounting Plate M35x0.5, also Latchable

- To insert unmounted optics with diameter 1/2", 1", 25 mm and 31.5 mm by means of retaining rings.
- Fitting retaining rings can be found in section Adapters.

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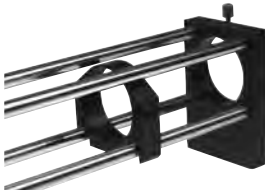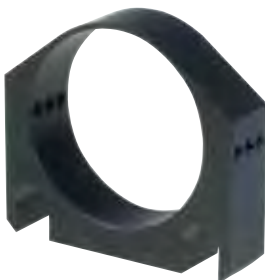Magnetic Cage Assembly Swapper Plates

The Magnetic cage assembly swapper plates allows the assembly of mounted optics into existing four-rod microbench cage assemblies. The swapper plate is equipped with two extra strong magnets made of neodymium-iron-boron (NdFeB). By that the plate is fixed at the desired position.

Also upside-down mounting is possible with a force up to 5N.


Two versions are available with diameter 25 mm and 30 mm.



Mounting plates with 10° angle to avoid back reflex.

- Strong magnets, magnetic force 5 N
- Can be used in four-rod assemblies
- For 25 mm and 30 mm mounted optics

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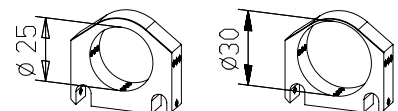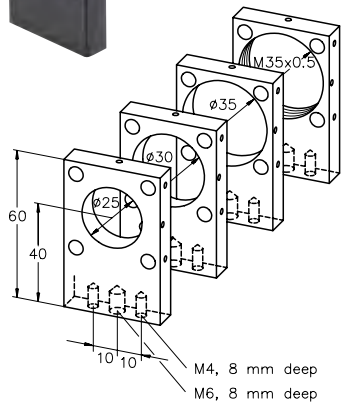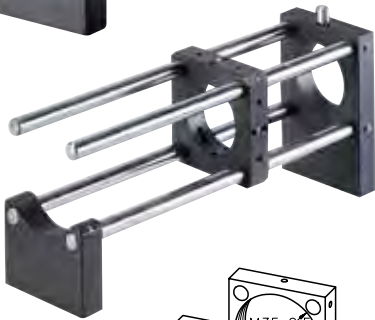
Holder 25/30/35/M35x0.5

- Holder for mounted optics with mount diameter of 25, 30 and 35 mm
- Holder M35x0.5 for unmounted optics with outer diameters of 1/2", 1", 25 mm and 31.5 mm with retaining rings
- Optical height of 40 mm
- Use with posts and holder possible
- Mounting options: fasten to mounting plates with M4 set screws, fasten to posts with M6 set screws, with base plate (e.g. G061673000) for freestanding/tabletop mounting
- Retaining rings for Holder M35x0.5 can be found in section "adapter" in this chapter.

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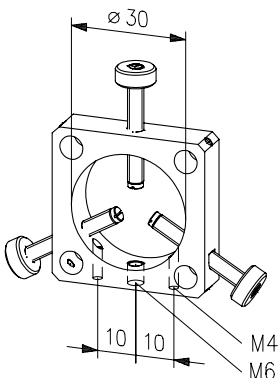Adjustable Holder 25

- For mounting cylindrical optical components from 3 to 25 mm diameter
- Mounting of components via screws
- Plastic tipped adjusting screws
- Gentle Clamping
- For longer components use two holders

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

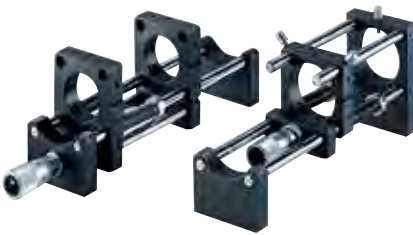
Rod Holder G/F/90°/V

The rod holder allows to mount microbench assemblies onto mounting plates, optical tables or posts.

- For assembling Microbench rods on breadboards
- Also mounts on posts
- Rod holder G: with a M10x0.5 tapped hole for accommodating a micrometer, i.e. G 10 (G061162000)
- Rod holder F: equipped with tapped-hole patterns, fasten from above with M6 or 1/4" screw
- Rod holder 90°: for right angle mounting of adapter plates,
- Rod holder V: through-hole M6, to be cut by yourself to the height you need for adapting to individual beam height
- Please see chapter positioning systems for micrometer screws

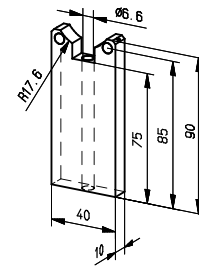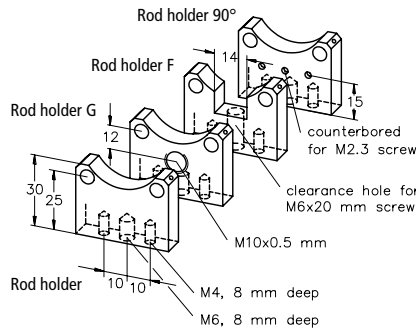Mounting example of Rod holder and Rod holder G with micrometer

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
Mounting Bases

These bases allow to mount microbench plates directly to an optical table or breadboard surface with M6 tapped holes on 25 mm centers or 1/4"-20 tapped holes on 1" centers. The optical axis height is 40 mm. The Mounting Base is fixed via a single screw and by this can be aligned at a desired angle.

- Direct mount of Microbench Plates on optical tables or breadboards with hole pattern 25 mm or 1"
- Mounting with screw M6 or 1/4"-20
- Mounting of the plate with screws M2.3x5
- Optical axis height 40 mm
- Delivery without Microbench Plate and screws

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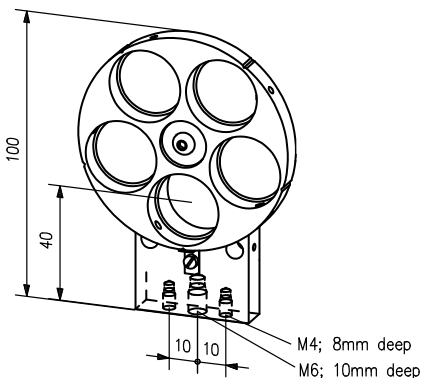Rotating Holder, 5 Positions

For mounting any construction element, in particular lenses, diaphragms and filters, with an outer diameter of 25 mm. The rotary disc is fixed in position by a detent spring. Two Rotating holders used in tandem create an ideal solution for use with neutral filters to achieve additional gradations.

- Accepts optics in \varnothing 25 mm mounts
- Spring-loading detents
- Various mounting options: mounts directly on \varnothing 6 mm rods, M4 tapped holes for mounting on bench plates, M6 central tapped hole for mounting on posts, mounts on tabletops with aid of mounting plates
- Threaded pins M2.3 included

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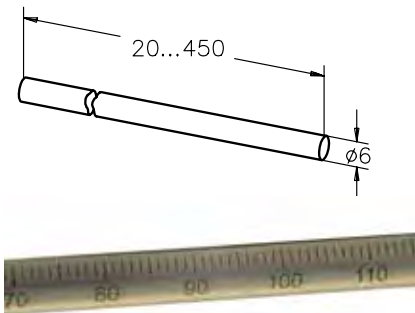
Rods


Rods are used to combine Microbench construction elements.

- Highest durability due to acid-resistant stainless steel
- Induction hardened -> no mark of screws etc. on the rods
- Lapped surfaces and tight tolerances guarantee optimum use with Microbench plates -> each part fits.
- Furthermore, the bending of these rods is absolutely minimum.
- Linearity 0.2 mm/m (linear decreasing)
- No tilting of the mounting plates.
- Rods 150 mm and 300 mm available with scale

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
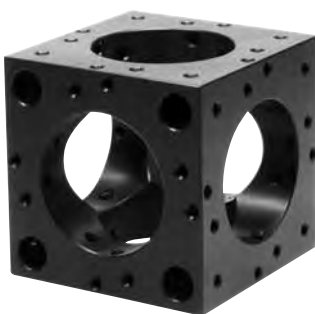
Cube

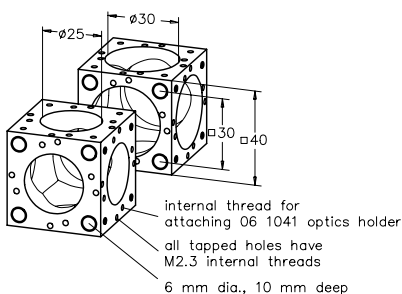- Multi-purpose structural elements
- For assembling compact, three-dimensional, optical systems, such as beamsplitters or beamsteering optics
- Cubes are delivered without screws, feel free to order: e.g. Set of screws M2.3 x 3 (qty. 150) G061011000
- Please look at section tools and screws regarding more accessories

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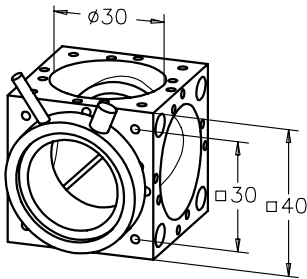Beam-Deflector Cube

Cubical mount equipped with rotatable plate holder

- Accepts 20x30 mm plane mirrors, beam-splitter plates, and right angle prisms
- Mirror reflecting surfaces aligned along axis of rotation
- Easy gluing by lateral fixation

- \varnothing 18 mm clear aperture
- Accepts G06 5092 000, positioning ring 30 (not included) for rotationally aligning optics


- Beam deflector cubes are delivered without screws, feel free to order: E.g. Set of screws M2.3x3 (qty. 150) G06 1011 000
- Please look at section tools and screws regarding more accessories

Umlenkwürfel Beam-Deflector Cube

Produktbezeichnung	Part No.
Item Title	Part No.
Beam-Deflector Cube	G061084000

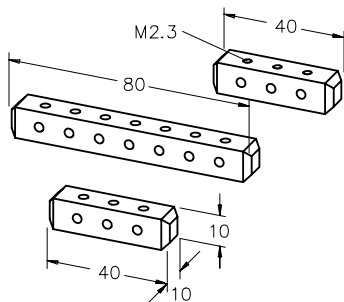Mounting example of Corner connector 40, in front you can see the Corner connector 40 with thread

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Corner Connector 40/80/40 with Threads

The Corner connector permits perpendicular connection of mounting plates to achieve orthogonal deflection of the optical axis or to split the beam.

- Corner connector 40 or 80 mm long for assembling mounting plates parallel or orthogonal to one another
- Corner connector 40 mm with threads for assembling components such as G061160000 Holder for micrometer

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
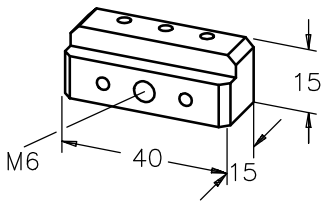
Corner Connector T

- For mounting plates on rods or breadboards
- Mounts on T 10 and TS 10 translation stages
- 40 mm length
- Please look at chapter "Positioning Systems" section "Linear Stages" regarding Linear stages T 10 and TS 10.

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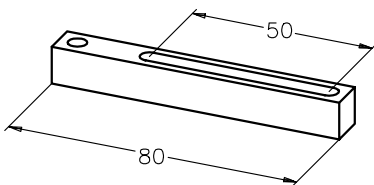Connector 80 V

The Connector 80 V for Microbench enables connections at various angles. It connects two mounting plates on a single plane. This Connector can be used to deflect the beam diagonally off axis. The slotted holes enable continuous adjustment of the angle and the mounting plates, and the crossed struts ensure optimum stability.

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- Links two coplanar mounting plates
- For adjusting distances and angles

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CAD files of all mechanical components can be found on the respective product sheets at www.qioptiq.com below the button "Docs and Drawings".

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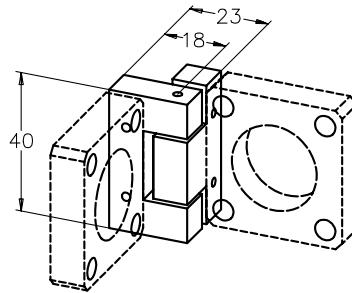
Swivel Connector

The Swivel Connector provides a backlash-free connection between two mounting plates, and has a pivot range of 120°. The ball-bearing mounted joint is pretensioned for backlash-free adjustment. The Connector 80 V is ideal for fixing the angle of the Swivel Connector.

- Combines pairs of mounting plates
- No-play precision swivel
- Pivots through 120°
- Use connector 80 V, G065068000, to fix swivel position


Gelenk

Swivel Connector

Produktbezeichnung	Part No.
Item Title	Part No.
Swivel Connector	G065067000

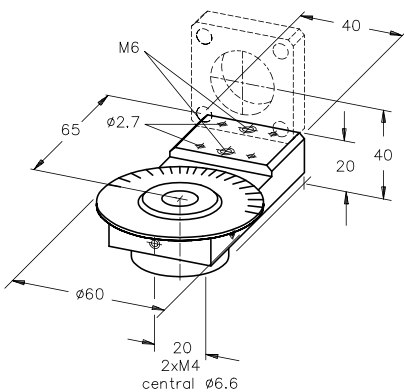Application example of Swivel arm with scale

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Swivel Arm with Scale

- For angular measurements in Microbench systems
- Accepts mounting plates or rod holder F
- Mounts on mounting plates, optical tables and carriers
- Optical axis height, 40 mm (when used with mounting plates)
- ± 90° scale with adjustable zero point
- Lockable with lateral clamping screw
- Mounting options: Central 6.6 mm Ø hole for top mounting, M4 threads (20 mm spacing) for bottom mounting

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Mounting example: Mounting bracket 40 at Post 25 with clamp

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


Mounting Bracket 40

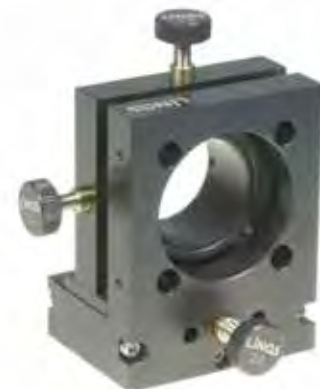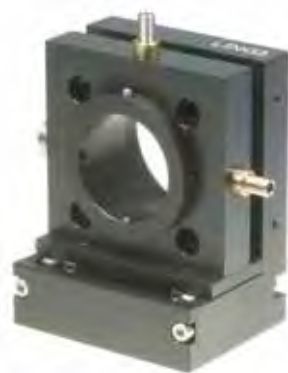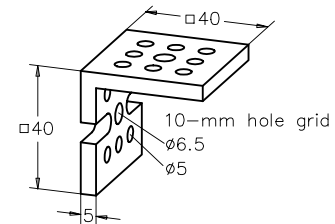- General purpose bracket for use in assembling Microbench compatible components
- Screws onto clamping fixture supplied with Post 25 with clamp
- Post 25 with clamp can be found in chapter "Mounts, Posts and Holders" section "Posts, Columns and Rods".

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XYZ-Fine Adjustment Unit

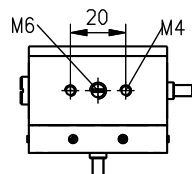
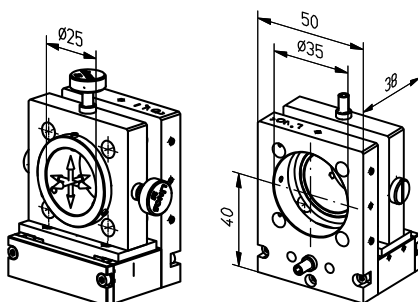
The XYZ-Fine adjustment unit enables a precise adjustment of components within the microbench cage system.

The inner diameter of 25 mm allows the use of Microbench mounted optics or elements like pinholes or fiber connectors.

The travel in XY-direction is ± 1 mm. In optical axis direction the travel range is 5 mm. Crossed-roller bearings ensure precise tiltfree guidance.

The XYZ-Fine adjustment unit is equipped with three stainless steel fine adjustment screws with 0.25 mm pitch. The screws have an internal hexagon socket, three knurled knobs will be delivered with the unit.

- Three fine adjustment screws with 0.25 mm pitch (100 TPI)
- Travel XY: ± 1 mm
- Travel Z: 5 mm
- Resolution: 1 μ m

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
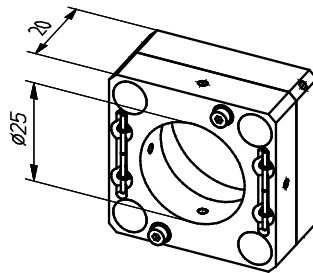
G06 1037 000

Slip Plate Positioner

The Slip Plate Positioner provides a coarse XY-adjustment of mounted optics in the microbench cages system.

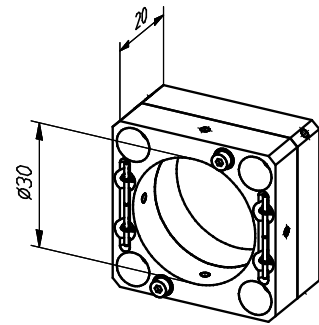- Two plates connected with springs
- Mounting diameter 25 mm and 30 mm
- Lockable
- ± 1 mm adjustment

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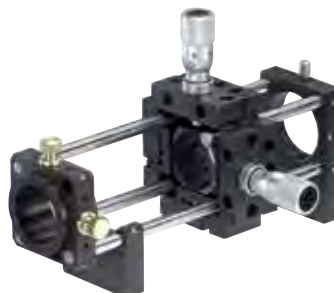
G061026000


Centering Mounting Plate/Holder 25

The Centering mounting plates and holders are compatible with any mount that has an outer diameter of 25 mm. Two fine-pitch screws are provided for adjustment of the centering fixture.

- Adjustable centering fixture
- High precision fine adjustment
- Inner ring rotatable
- ± 1 mm travel
- Screws with 0.25 mm pitch
- Threaded screws M2.3 included

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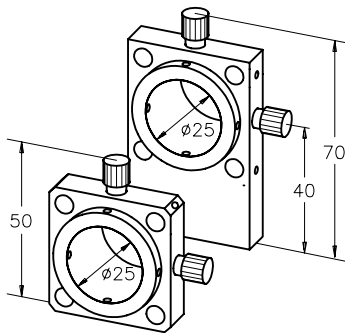
X-Y Precision Translation Stage

The XY-adjustment is available in two versions, with micrometer screw or with fine adjustment screw.

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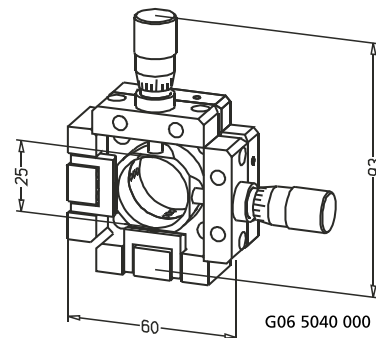
- Accepts Ø 25 mm components
- Micrometer screws and fine-adjustment screws with 0.25 mm pitch
- 0.01 mm cylinder scale
- ±1 mm adjustment range
- Connection thread of the screws M10x0.5
- Threaded pins M2.3 included
- G06 5070 000 with fine adjustment screws, otherwise identical to G06 5040 000

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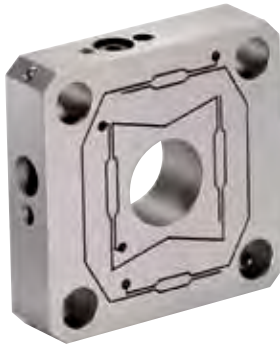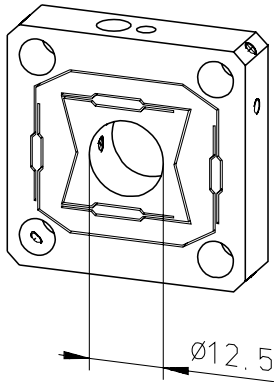
fle.X-Plate

The fle.X-plate is a monolithic adjustment unit with flexures.

It is available in different versions:

- fle.X-plate XY Al and fle.X-plate XY Steel permit an adjustment range of ± 0.25 mm. Adjustment is carried out using tension and pressure screws, which are subsequently secured in position with lock-nuts for excellent long-term stability.

- The fle.X-plate XY Diff has differential screws for high-precision adjustment on two axes. The differential screws have knurled knobs for easy adjustment. Material: Aluminum black anodized.


The pitch is 0.1 mm per rotation.

- For fiber optics, spatial filters, pinholes
- Flexure from hardened steel or aluminum
- Excellent long-term stability

- Accepts components with outer diameter 12.5 mm
- fle.X-plate XY Al/Steel, spindle pitch: 0.50 mm
- fle.X-plate XY Al/Steel, reproducibility: very good, since play is minimized by tension and pressure screws
- fle.X plates are delivered with tools for adjustment screws.
- For fixing in the Microbench you will need screws M2.3, you can order the following: Set of screws M2.3x3 (qty. 150) G06 1011 000.

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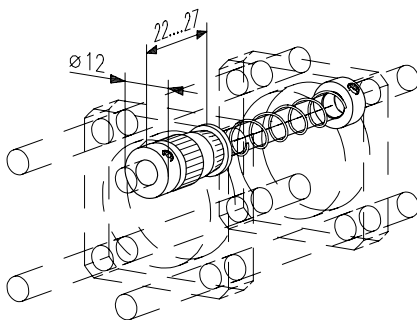Z-Adjustment M

The Z-Adjustment M, consisting of fine-adjustment screw, compression spring and collar is a highly precise fine-adjustment with 5 mm translation range. It allows a backlash-free adjustment of optics, pinholes, fiber plugs etc., has a very compact design and later integration into existing setups is possible.

- High precision fine-adjustment with 5 mm travel
- Consisting of fine-adjustment screw, compression spring and collar
- Backlash-free adjustment of optics, pinholes, fiber plugs etc.
- Compact
- Later integration possible

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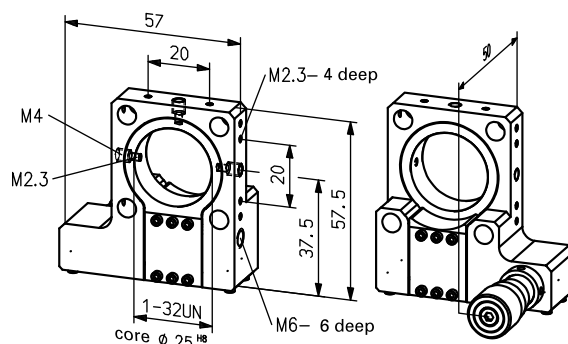
Z-Fine Adjustment M

This is the most precise Z-fine adjustment of our Microbench elements with a direct mount for 25 mm outer diameter components. The Z-Fine Adjustment is equipped with a fine-adjustment screw with 0.25 mm pitch. The translation range is 1.5 mm with an accuracy of 1 µm. Because of the 1-32UN internal thread this unit fits to our Tube Mounting System C. This component is ideally suitable for example for fiber alignments due to its exact centering without nearly no tilt.

The Z-Fine Adjustment is available in two versions, with micrometer or with fine adjustment screw. The micrometer screw has a pitch of 0.25 mm. The spindle has a hexagon socket and can be driven by the provided knurled hexagon knob. If the hexagon socket is not needed, the knob can be glued in.

- High precision fine-adjustment
- Clearance-free adjustment of optics, pinholes, fiber plugs, etc.
- For components with outer diameter of Ø 25 mm
- Compact
- 1.5 mm travel
- Micrometer screws with 0.25 mm spindle pitch
- Accuracy 1 µm
- Connection thread of the micrometer M10x0.5
- With 1-32UN inner thread (C-Mount)
- Micrometers are described in chapter "Positioning Systems"

Z-Feinverstellung M Z-Fine Adjustment M

Produktbezeichnung	Part No.
Item Title	Part No.
Z-Fine Adjustment M with Micrometer	G061061000
Z-Fine Adjustment M with Fine Adjustment Screw	G061063000

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Linear Stages for Microbench

The linear stages are available in various versions, central or offset drives, with micrometer screw or with fine adjustment screw.

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- Translation stages for alignment along Z-axis of components in mounting plates for assembly into Microbench
- Choice of central (type T) or offset drives (type TS)

- 10 mm travel
- 10 µm positioning resolution
- Mounting of stages via corner connector T, G06113000 (2 pcs. per stage)

- Adaption to Microbench mounting plates with the help of corner connector T, require 2 ea. per translation stage for installation.
- Please look at section Positioning Systems, Linear stages regarding technical data and drawings.

T 10 M

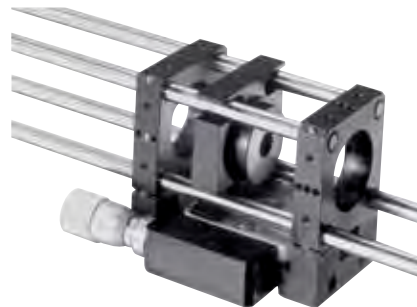
TS 10 M

T 10 F

TS 10 F

Mounting example of linear stage TS 10 M with 2 corner connectors T

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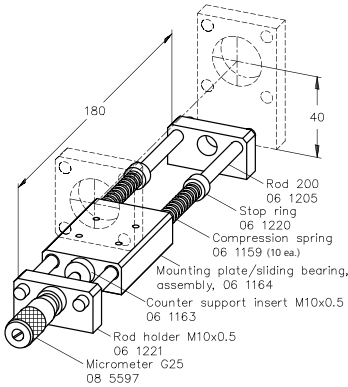
Z-Translation Stage with Micrometer

- Manual translation stage for optical axis height of 40 mm
- Direct mounting of Microbench mounting plates
- For assembling on mounting plates or to integrate into rod structures
- All components are available separately

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- With micrometer, readable to 0.01 mm
- Connection thread of the micrometer M10x0.5
- 25 mm travel

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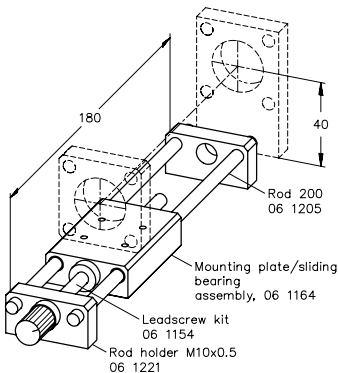
Z-Translation Stage with Fine-Adjustment Screw

- Manual translation stage for optical axis height of 40 mm
- Direct mounting of Microbench mounting plates
- For assembling on mounting plates or to integrate into rod structures
- All components are available separately

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- Fine-adjustment with 0.5 mm screw pitch
- 105 mm travel

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
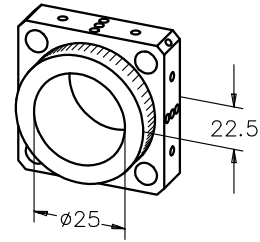
Rotary Support

- For mounting slits, reticles, polarizers, prisms etc., requiring rotational alignment
- Angular scale graduated in 5° increments
- Rotation axis lockable with Screw M2.3

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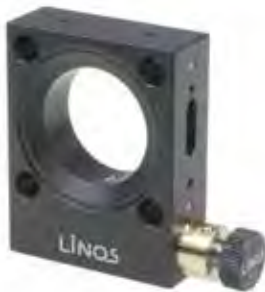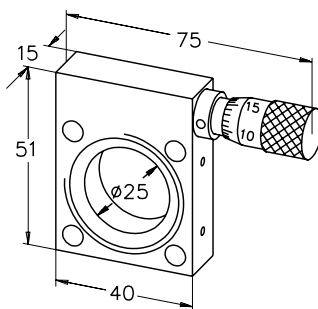
Rotary Support 20°

The rotary support is available in two versions, with micrometer screw or with fine adjustment screw.

The fine-adjustment screw with 0.25 mm pitch has a hexagon socket and can be driven by the provided knurled hexagon knob. If the hexagon socket is not needed, the knob can be glued in.

- Ideal for mounting polarizers and birefringent filters in laser technology
- Micrometer version equipped with Micrometer Screw G10
- Angular-adjustment range ±10°
- Angular sensitivity 10"
- Connection thread of the micrometer M10x0.5
- Ø 25 mm clear aperture

Drehfassung 20° Rotary Support 20°

Produktbezeichnung	Part No.
Item Title	Part No.
Rotary Support 20° with Micrometer Screw	G065088000
Rotary Support 20° with Fine-Adjustment Screw	G061064000

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The Motorized Rotation Mount enables fast rotation of components like polarizers. There are two versions, with servo motor or stepper motor available.

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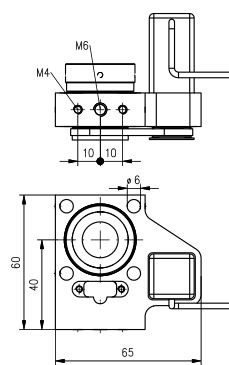
The measurement of the rotation speed is possible with the optional photo sensor.

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
- *For fast motorized rotation, periodic modulation*
- *Servo- or stepper motor*
- *Positioning of optical components (stepper)*
- *Optional photo sensor*

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- *For elements with 25 mm diameter*
- *Up to 400 rpm*

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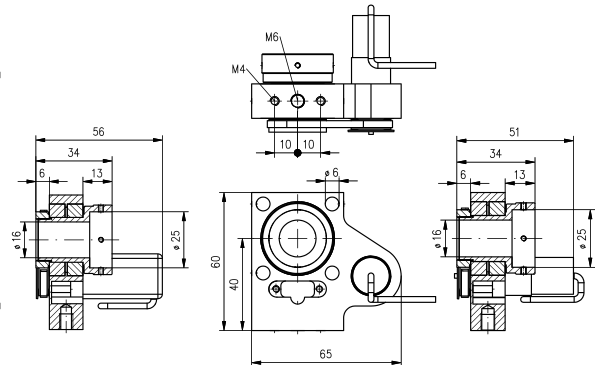
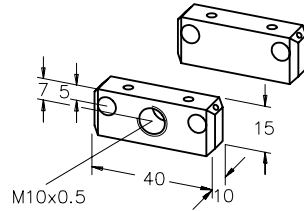
In front: Holder for Micrometer

Counter Support

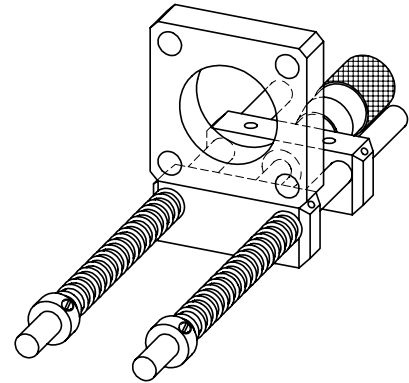- For use with G06 1160 000 holder for micrometer
- To realize a Z-translation

Example for a Z-translation with Holder for Micrometer, Micrometer G 10, Counter support, Compression springs and stop rings

Gegenlager Counter Support

Produktbezeichnung	Part No.
Item Title	Part No.
Counter Support	G061161000

Example for a Z-translation with Holder for Micrometer, Micrometer G 10, Counter support, Compression springs and stop rings

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

Holder for Micrometer

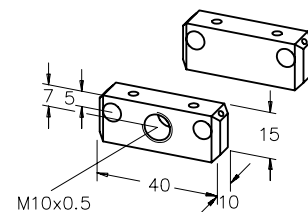- Fits to mounting plates
- Mounts on Microbench rods
- Ideal to realize a Z-translation with Micrometer G 10, counter support, compression springs and collars
- For micrometer with thread M10x0.5
- Please see chapter Positioning Systems, section "Micrometers and Adjustment screws" for micrometer screws

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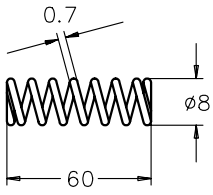Compression Spring, 60x8x0.7

- Ideal to realize a Z-translation with Micrometer G 10, counter support, compression springs and collars
- Slips over \varnothing 6 mm rods
- Two Compression springs are needed for assembly of a Z-translation stage.

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Adjustable Prism Insert 30

- Surface for cementing prisms or beam splitter cubes with edge length of 20 mm
- High precision fine adjustment
- Three screws for adjusting tilt angles and setting axial position
- Fits into G06 1081 000 Cube 30
- Rotatable in mounting plate

- Fine-adjustment screws with 0.15 mm pitch
- Screws with internal hexagon
- Two Allen wrenches included
- Independently adjustable axes
- \varnothing 8 mm (clear aperture) through bore

- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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Please find detailed information about glueing optics on prism and mirror holder in our Brochure "Microbench - Introduction, solutions and applications" (download at www.gioptiq.com). You will find the brochure for download at each Microbench product under "Docs + Drawings"

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
Adjustable Mirror Insert 30

- Surface for cementing mirrors or beam splitting plates with size 22.4 x 31.5 mm
- High precision fine adjustment
- Rotatable in mounting plate
- Fits into G06 1081 000 Cube 30
- Fine-adjustment screws with 0.15 mm pitch
- Screws with internal hexagon
- Two Allen wrenches included
- Independently adjustable axes
- \varnothing 14 mm (clear aperture) through bore in mirror mounting surface
- \varnothing 8 mm (clear aperture) through bore the base
- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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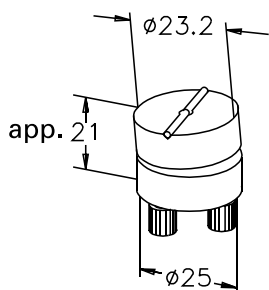
Tilting Platform

- Flat surface for cementing prisms, mirrors, beam splitting cubes, etc.
- Fits between pairs of Microbench rods
- Two thumbscrews adjusting angular alignment by $\pm 4^\circ$
- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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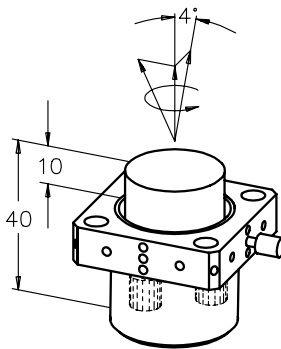Prism Table 25, Adjustable

- Prism mount is rotatable in mounting plate
- High precision fine adjustment
- 3 screws for adjusting tilt angles and setting axial position
- Components cement into place
- Protective cap for preserving thumbscrew settings

- High precision fine adjustment (screws with 0.25 mm pitch)
- Three thumbscrews adjusting angular alignment by 4° in all directions

- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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Mounting Tool for Beamsplitter Cubes

The mounting tool will help you to mount beamsplitter cubes or prisms centric on the adjustable prism insert 30 (G063731000) or prism support G065089000. The tool is made from plastic not to damage the surface of the optics.

- Two sizes, for beamsplitter cubes 16 mm or 20 mm
- Also useable for 90° prisms
- Material black plastic

- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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Rotation Adjustment for Cube Inserts

The rotation adjustment for cube inserts allows the precise adjustment of prism or mirror inserts with diameter 30 mm in microbench cubes 30.

The adjustment is equipped with a fine adjustment screw with 170 TPI and allows an adjustment range of $\pm 2.5^\circ$.


The screws have an internal hexagon socket, a knurled knob will be delivered with the unit. If the hexagon socket is not needed, the knob can be glued in.

The adjustment unit will be mounted on one side of a cube 30. All cube inserts with diameter 30 can be used.

- Precise angle adjustment of cube inserts
- For Cube inserts 30 mm diameter
- Easy mounting
- Fine-adjustment screws with 0.15 mm pitch
- Screws with internal hexagon

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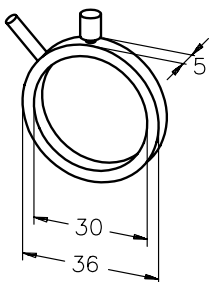
Positioning Ring 30

- Rotational alignment of $\varnothing 30$ mm mirrors, prism mounts etc.
- Equipped with adjusting lever

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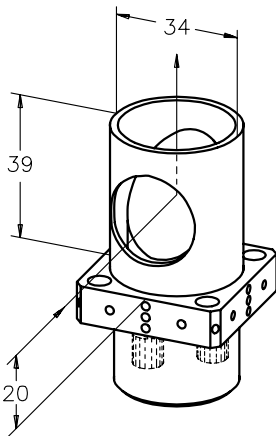Beam Steering Mirror Assembly


- Elliptical (22.4 x 31.5 mm) front surface plane mirror (G340003000) inclined at 45° in mounting plate
- UV type with UV plane mirror (G340094000)
- High precision fine adjustment
- Three screws for adjusting alignment and setting axial position
- Rotatable through 360°
- Protective cap for preserving thumbscrew settings

- Screws with 0.25 mm screw pitch
- $\pm 4^\circ$ alignment range

- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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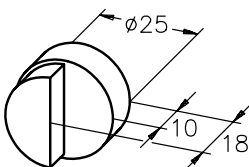Optic Support 25



Diameter 25 mm. The optic support 25 allows to use unmounted mirrors, beamsplitter plates, filters etc. into the beam path. The elements will be glued against the grade, that the surface will be in the center of rotation.

- For inserting unmounted mirrors, beamsplitter plates, filters, etc. into beam path
- Optical components cement into place

- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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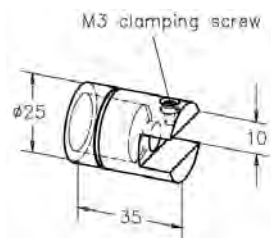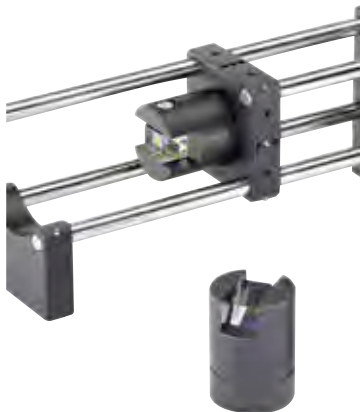Clamp Holder for Prism

- For mounting optical components (prisms, beam splitting cubes, thin-film polarizers, 10x17x 8.5 mm neutral density wedge filters, etc.) with 10-mm edge lengths
- Polyamide screw provides gentle clamping action
- Central bore allows inserting mount in beam path

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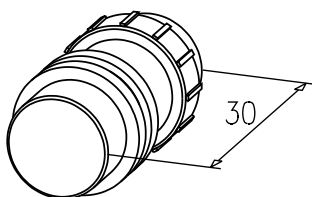
Prism Support 25


The Prism support consists of a threaded bolt with a 25 mm retainer ring and locknut, and is used to mount deflector prisms, beam-splitter cubes, 30° prisms or similar components in the beam path. The optical construction elements are either bonded onto the end plane of the threaded bolt or clamped between two threaded bolts.

- For mounting beam splitting cubes, prisms, etc.
- Optical components may be clamped between two supports or cemented into place
- Threaded bolt with a 25 mm retainer ring and locknut.
- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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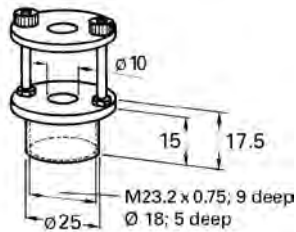
Prism Table 25

- For mounting prisms, plane plates, or beam splitting cubes
- With central bore for beam passage
- 20 x 30 mm unobstructed mounting surface
- Max. component height accommodated is roughly 25 mm

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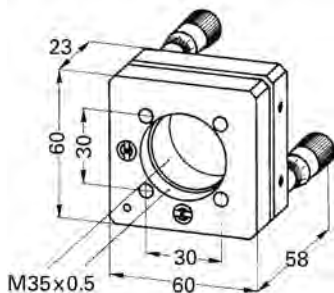Adjustable Optic Holder 31.5

The Adjustable Optic Holder is available in two versions, with micrometer screws or with fine adjustment screws.

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- Precision, high-resolution cinematic optic mount
- With micrometers (type M) or ungraduated fine-adjustment screws (type S)
- Includes two retaining rings (G065069000) for diameter 31.5 mm optics
- Accepts Ø 25, 1", 1/2" optics with suitable retaining rings
- ± 4° angular alignment range
- Micrometer drives with 0.8' angular-resolution
- 0.25 mm lead screw pitch
- Accommodates optical components with a thickness of 1.5 to 10 mm
- M35x0.5 internal threads on both plate apertures
- Mounting options: M6 tapped hole in base for mounting on posts or mounting plates, mounts directly on Microbench rods
- See the "Adapters" section for details on threaded rings for unmounted optics with an outer diameter of 25 mm, 1" or 1/2".

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
Overview Table Adapter

This is an overview on the adapters of Microbench, Nanobech and Tube system. Detailed information on these adapters can be found on the following pages and in the respective chapters.

Außen-Ø (mm)	Innen-Ø (mm)	Länge (mm)	Artikel-Nr.
<i>Outer ø (mm) respectively thread</i>	<i>Inner connector respectively ø (mm)</i>	<i>Length (mm)</i>	<i>Order-No</i>
12.5	FC	8	G06 7023 000
12.5	SMA	8	G06 7024 000
16	FC	8	G05 0310 000
16	SMA	8	G05 0311 000
25	23.2	10	G03 8855 000
25	16	10	G05 0305 000
25	15	12	G06 1626 000
25	25	20	G06 5042 000
25	FC	10	G06 7054 000
25	SMA	10	G06 7055 000
30	C-Mount	10	G03 8851 000
30	25	16	G06 1602 000
30	23.2	20	G06 1634 000
30	C-Mount	16	G06 1658 000
30/1.035-40UN	M23.2x0.75	13	G06 0407 000
30	1.035-40UN	10	G06 1168 000
35	30	16	G06 1637 000
35	30, 3.5° tilt	16	G06 1639 000
50	For Mounting plate	13	G06 1674 000
25/30	Full material	20	G06 1601 000
25/30	W0.8"x1/36"	20	G06 1621 000
25/30	16	20	G06 1622 000
25/30	19	20	G06 1623 000
25/30	23.2	20	G06 1636 000
25/30	C-Mount	22	G06 1664 000
25/C-Mount	16/W0.8"x1/36"	34	G06 1661 000
30/C-Mount	M19.5x0.5	34	G06 1669 000
C-Mount	Connection for Microbench rods	21	G06 1659 000
C-Mount	M23.2x0.75	8	G06 1667 000
C-Mount	W0.8"x1/36"	8	G06 1668 000
M17x0.75	C-Mount	11	G03 8878 000
M35x0.5	30	16	G06 1638 000
W0.8"x1/36"	C-Mount	14	G06 1662 000
30 (e.g. for Cube 30)	SMA Collimator		G06 1087 000
Connection for Microbench rods	SMA Collimator		G06 1088 000
25 (e.g. for Cube 25)	Full material	5	G06 1079 000
30 (e.g. for Cube 30)	Full material	5	G06 1078 000
40x40, Adapter plate for FLS 40 carrier	Full material		G06 1085 000

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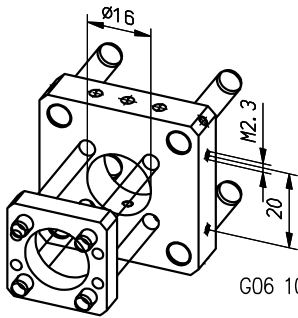
Microbench - Nanobench Adapter

To connect the cage mounting systems Microbench and Nanobench.

- Rod holes for Nanobench rods, hole pattern 16 mm
- Rod holes for Microbench rods, hole pattern 30 mm
- Inner diameter 16 mm for optics mounted in Nanobench mount
- The plate will be delivered without screws. Please use M2.3 Microbench and M2.5 Nanobench screws to fix the rods.

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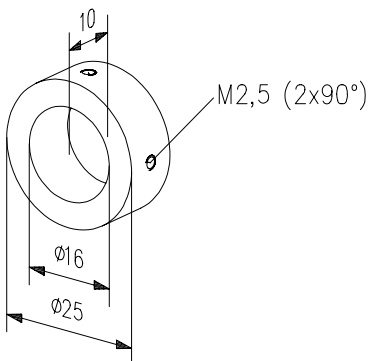
Adapter 16/25

To fit mounted Nanobench optics (outer diameter of 16 mm) in Microbench components with internal diameter of 25 mm, e.g. Mounting Plate 25, G06 1010 000

- To fix the mounted lenses, you will need Set of Setscrews M2.5x3, (G05 1520 000), not included

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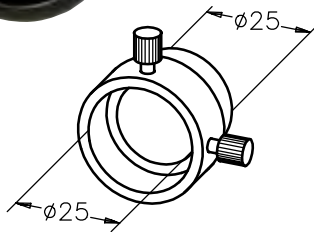
Adapter Collar

Allows rapidly interchanging of components with mounts outer ϕ 25 mm or tube ϕ 25 mm

- Two M2.3 thumb screws for mounting components are included

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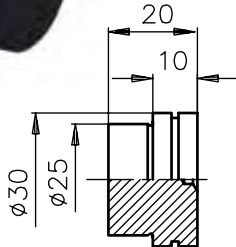
Cap 30/25

- For closing off ϕ 25 or 30 mm apertures at ends of beam paths
- For fabricating custom optic holders and other special components

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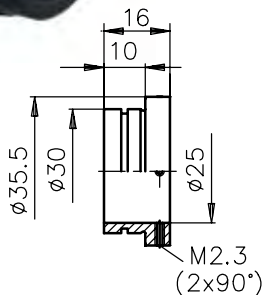
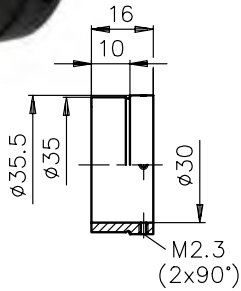
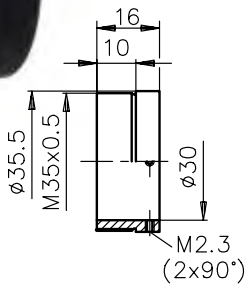
Adapter M35x0.5/30

For mounting Ø 30 mm components on items with M35x0.5 threads

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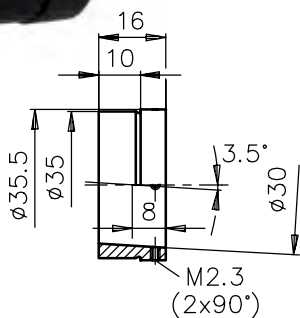
Adapter 35/30-3.5°

• Optical components are fixed at an angle of 3.5° to avoid backreflections

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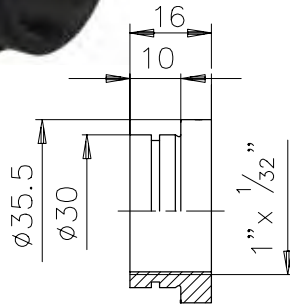
Adapter 30/C-Mount

- For mounting devices, such as CCD cameras, laser beam expanders, etc., with C-Mount (1" x 1/32) threads in Ø 30 mm apertures

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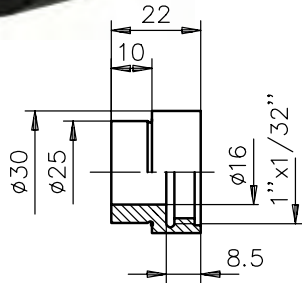
Adapter 25/C-Mount

- Outside diameter 25 mm and internal C-Mount thread

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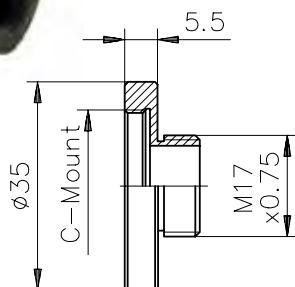
C-Mount-Adapter-M17

- For fitting objectives to C-Mount threads

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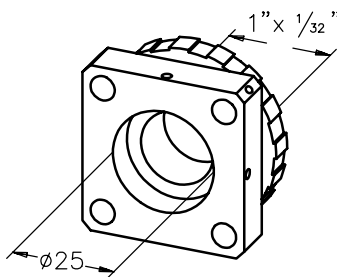
Adapter C-Mount/Microbench

- For mounting lasers or CCD cameras with C-Mounts to Microbench assemblies
- Knurled retaining ring for locking alignment at the C-mount thread
- Fits directly on Microbench rods
- Diameter 25 mm central aperture

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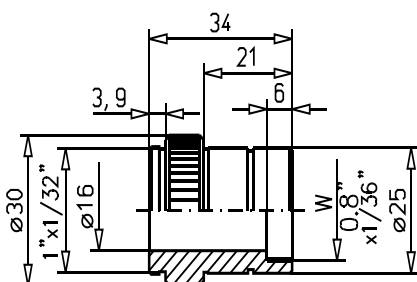Adapter 1" to W0.8" or Ø 25 mm

- For mounting components with C-Mounts on standard microscope threads
- Fits to mounting plates

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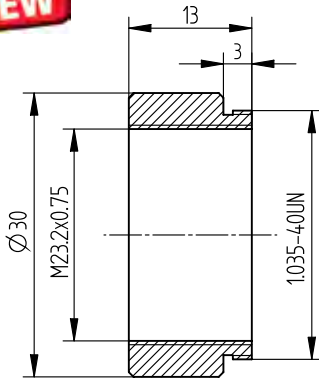
Adapter Bench Systems 30 & 1.035''-40 UN


For adapting other bench/ cage systems with the LINOS Microbench. One side fits to Microbench mounting plates with 30 mm diameter, the other fits into mounting plates with 1.035''-40 UN internal threading. Furthermore, the adapter has an M 23.2 x 0.75 internal threading.

- Combines various cage systems
- Adapts mounts and carriers designed for imperial units to Microbench standard
- Suitable for Microbench Mounting Plates Ø30 and plates with 1.035-40 threading
- Internal thread allows mounting of optics up to Ø22.4 via retaining rings

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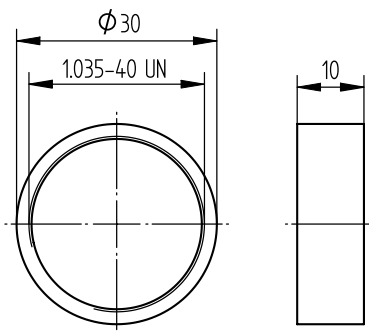Adapter 1.035''-40UN / 30

For adapting mounted optics with 1.035''-40 UN threading to Microbench mounting plate with 30 mm diameter.

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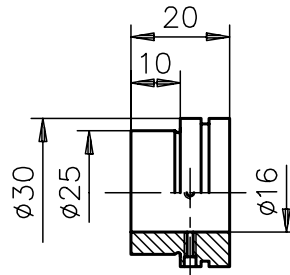
Adapter 16

- Fits outer- \varnothing d = 16 mm compact objectives to \varnothing 25 and 30 mm mounts

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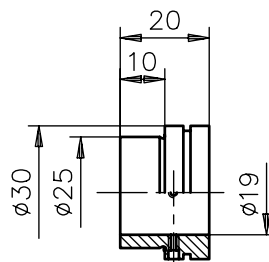
Adapter 19

- Fits outer- \varnothing d = 19 mm compact objectives to \varnothing 25 and 30 mm mounts

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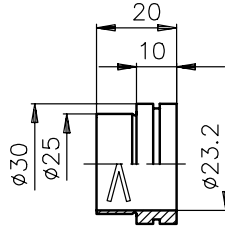
Adapter 23.2

- For mounting microscope eyepieces with standard Ø 23.2 mm in mounting plates 25 or 30

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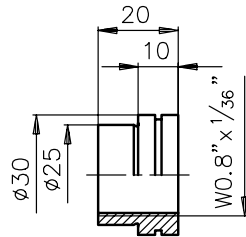
Adapter W0.8''x1/36

- For mounting microscope objectives with standard W0.8''x1/36" (RMS) threads

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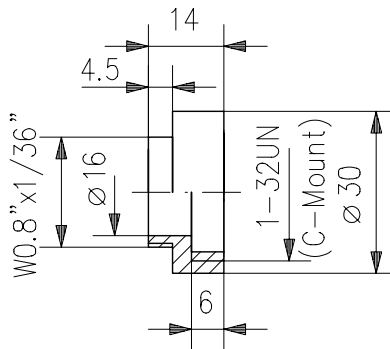
Adapter W0.8'' to 1''

- For mounting components with C-Mount threads onto standard microscope threads.

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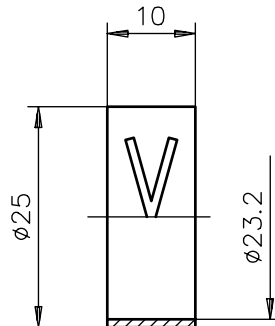
Adapter 23.2 to Ø 25

- For mounting microscope eyepieces with standard Ø 23.2 mm in Ø 25 mm apertures

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
Produktbezeichnung	Part No.
Item Title	Part No.
Adapter 23.2 to Ø 25	G038855000

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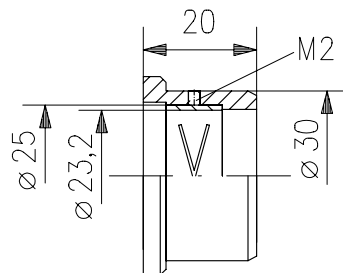
Eyepiece Adapter 23.2/30

- For mounting microscope eyepieces with standard Ø 23.2 mm in Ø 30 mm apertures
- Tube length retains unchanged after adaption

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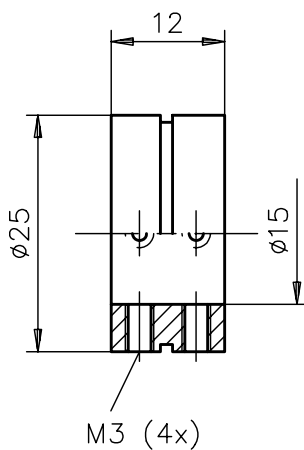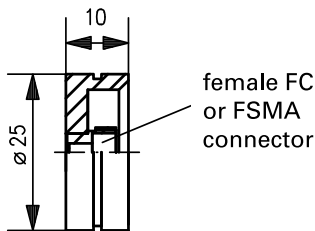
FC/FSMA Fiber Adapter

- *Mounting flanges without optics*
- *Accept fibers with FC connector / FSMA connector*
- *Outer Ø of 25 mm fits to mounting plate 25, holder 25, centering mounting plate 25 etc.*

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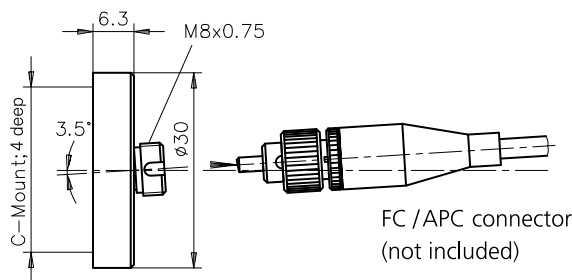
FC Fiber Adapter 3.5°

- *Mounting flange without lens*
- *For use with FC / APC and AFC connectors*
- *Compensates for beam deflection that occurs when FC / APC and AFC connectors are used*
- *C-Mount female thread*
- *Outer diameter of 30 mm for assembly on Microbench Mounting plate 30 or Holder 30*

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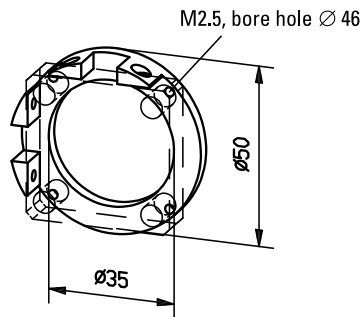
Instrument Flange

- Adapts microbench mounting plates to various instruments housings

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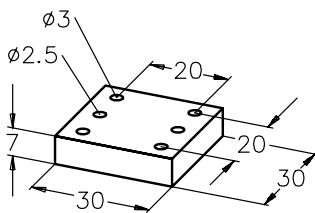
The instrument flange (G06 1674 000) enables connecting external modules to your system, e.g. for laser beam coupling in a microscope. View further information in our brochure "Microbench, Bench and Rail systems - Introduction, solutions and applications". You will find the brochure for download at each Microbench product under "Docs + Drawings".

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

Adapter Plate

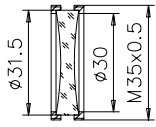- For mounting Microbench plates on e.g. a TB 50-16 positioner

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
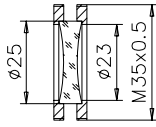
Set of Retaining Rings

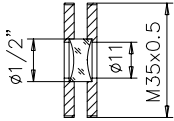- Set each 2 retaining rings
- To mount optics diameter 31.5 mm, 25 mm, 1/2" or 1" in mounting plate M35x0.5 or holder M35x0.5

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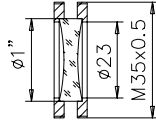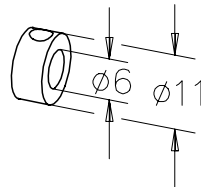
Set of Collars

- For use as clamps or adjustable stops on Microbench rods
- To realize a Z-translation in combination with rods, counter support, holder for micrometer, micrometer G 10, and compression springs

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Two Beam Interferometer

Many physical occurrences can be described in terms of waves. The emission of radio waves and the emission of light waves from light sources are all examples of this phenomena. If we overlap these waves, we can observe the phenomena known as interference. It is sometimes much easier to view the desired interference effects by introducing a disturbance into the process. Often, we carefully guide the interference to obtain highly accurate measurements of physical sizes. These physical sizes are determined by the path difference of numerous light beams (two in this case). Distances, angles, velocity, geometrical forms, flatness and many other measurements can be determined by interference and compared to the wavelength of the light source.


In turn, an interferometer with known geometrical data can be configured to determine the wavelength of a light source and analyze the spectral characteristics of this source.

This interferometer, assembled out of Microbench components, can be used to study the basic principles of two beam interference, perform exact measurements and can be expanded for other applications.

- Rigid setup
- To be combined with laserset (G064016000, please look at section microbench sets). Suitable laser are described in chapter "Instruments", section "Light Sources and Laser".

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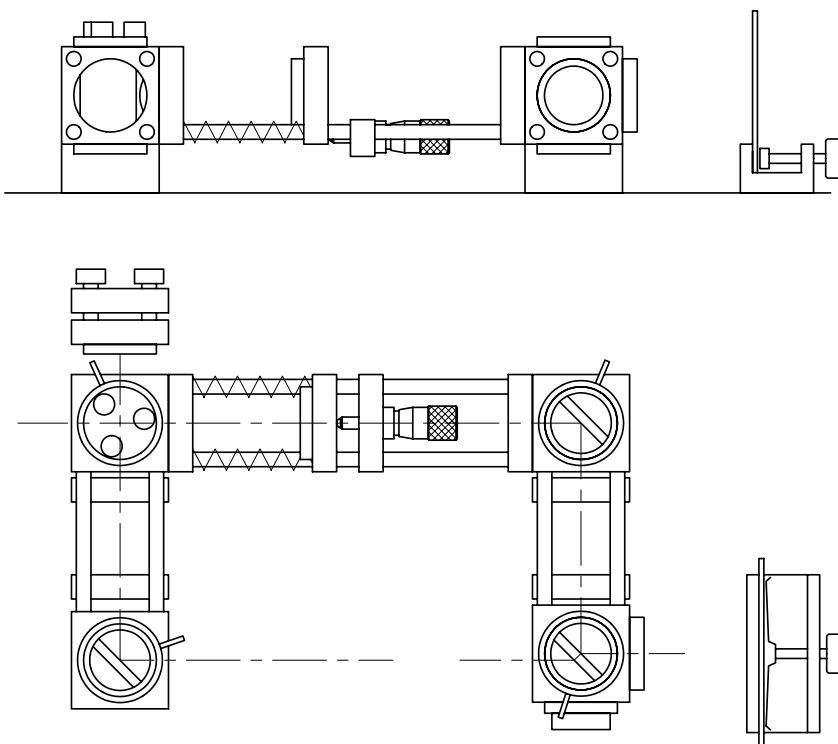
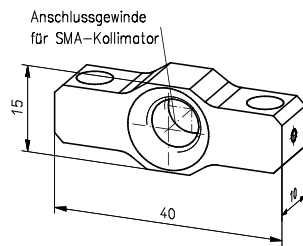
Collimator Holder

Collimators with SMA thread can be integrated into the microbench system with the collimator holder.

The Collimator holder S has two holes for the microbench rods so it can be positioned at any place on the rods.

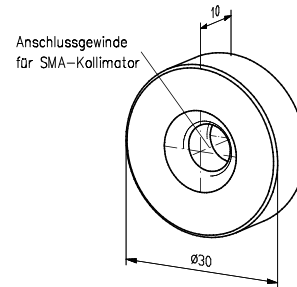The Collimator holder W has an outer diameter of 30 mm to be integrated into microbench cube or mounting plate 30.

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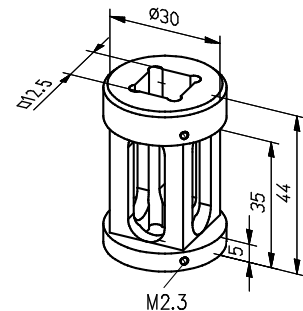
Cuvette Holder

The Cuvette holder can be integrated into the cube 30 and is used to hold standart cuvettes with edge lengths of 12.5 mm.

The cuvette holder will be delivered with one cover cap.

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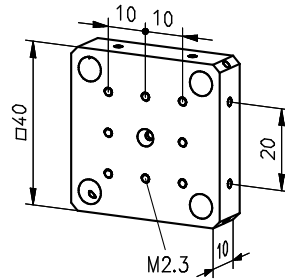
Mounting Plate FLS

The Mounting plate FLS can be used for vertical mounting of microbench setups at carriers of the FLS 40 system.

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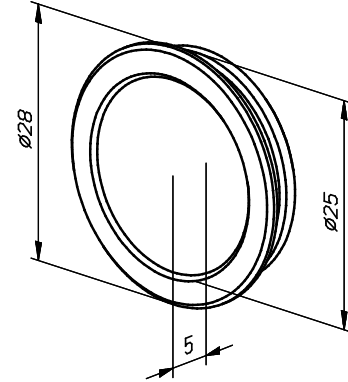
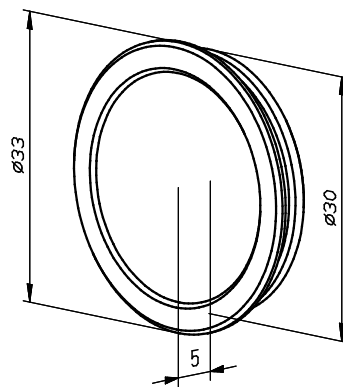
Closing Caps

The Closing Caps can be used to close unused apertures in microbench cubes. There are two sizes with diameter 25 and 30 mm available. The caps are made of black anodized aluminum.

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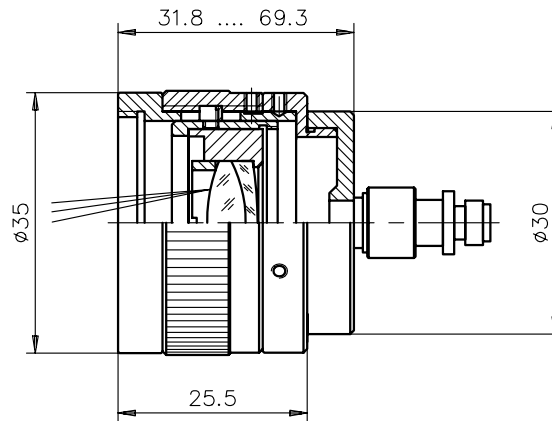
Focusing Collimators

- To focus or collimate divergent light
- Compact design
- Excellent beam quality
- Low wavefront distortion
- Achromats with antireflection-broadband coating
- 5 mm travel
- For spectral range 400 - 800 nm wavelength
- For singlemode fibers with NA = 0.1 or multimode fibers with NA = 0.22
- Two FC connector versions: for fibers with 8° polish (FC8/PC8) or for fibers with 0° polish (FC0/PC0)

Focusing collimator with achromatic lens

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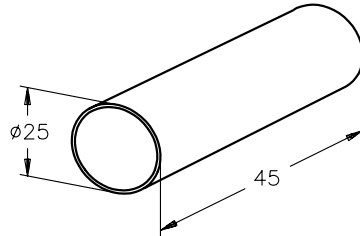
Tube

- Blocks ambient / scattered light
- May also be used to hold eyepieces
- Outer-Ø 25 mm, inner-Ø 23.2 mm

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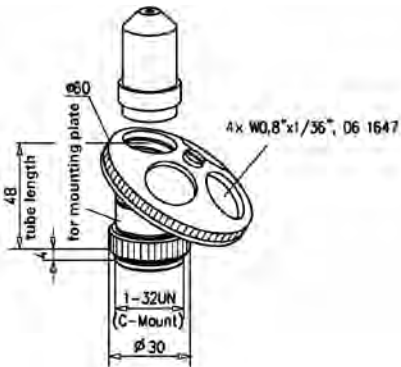
Lens Turret 4x with C-Mount Adapter

- For mounting 4 microscope objectives
- Spring-loaded detents
- Removable C-Mount adapter
- To combine with Mounting plate 25
- Ball-bearing turret disk Ø 60 mm
- Fitting threads:
 - With adapter: 1-32 UN (C-Mount)
 - Without adapter: W0.8"x1/36"

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Microbench Alignment Plate

The Microbench Alignment Plate will help you to adjust the beam path into microbench assemblies. The plate is made from anodized aluminum, with a hole diameter 1 mm at the optical axis and printed crosshairs.

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Circlip Kit

- Set of circlips mating to grooved reduction rings
- Consists of 5 circlips for Ø 25 mm reduction rings, 25 circlips for Ø 30 mm reduction rings plus a pair of circlip pliers

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Special Screwdriver

- For threaded pins M2.3
- To be used in small space mountings
- 1.8 mm blade width
- 20 mm shaft length

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Screwdriver

- To screw threaded pins M2.3 in mounting plates, holder, rod holder or cubes
- 1.8 mm blade width

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Set of Threaded Pins M2.3x3

- Accessories to mounting plates, holdes, cubes etc.
- To clamp components onto rods
- Set = 150 pieces

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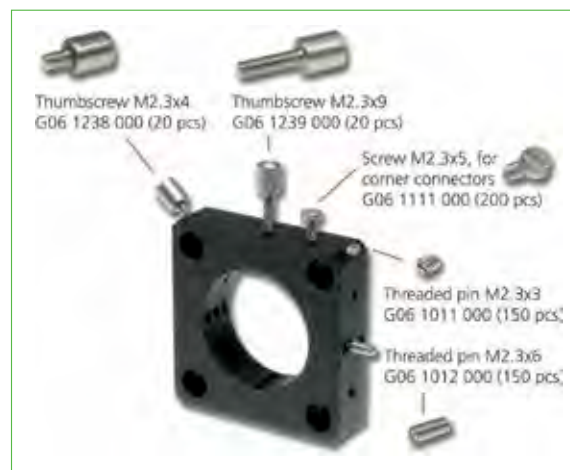
Set of Threaded Pins M2.3x6

- To fix mounted optics in mounting plates and holders
- Set = 150 pieces

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Set of Thumbscrews M2.3x9

- To fix mounted optics in mounting plates 25/30 and holders 25/30
- Set = 20 pieces

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Set of Thumbscrews M2.3x4

- To clamp mounted components into mounting plates 35 or holder 35
- For quick fixing of mounting plates, holders and rod holders on rods
- Set = 20 pieces

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Set of Screws M2.3x5


- To mount corner connectors to mounting plates
- Set = 200 pieces

Satz Schrauben M2.3x5 für Eckverbinder Set of Screws M2.3x5 for Corner Connector

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Set of Screws M2.3x5, (200pcs.)	G061111000

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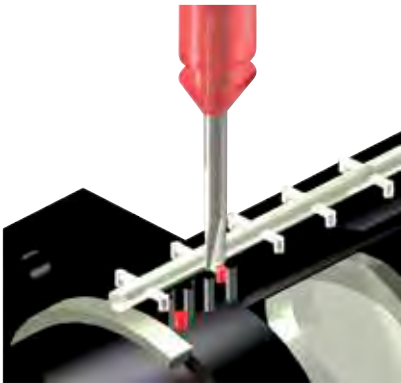Polyamide Plugs

- *The plugs prevent optics from tipping over during rotation or subsequent attachment (no backlash)*
- *The polyamide plugs are inserted into the thread before the setscrews M2.3 are attached*
- *1.5 x 1.5 mm (qty. 100, 10x10)*

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*Mounting and production of mechanical
components in Goettingen about 1910*

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*Setting of parameters at one of our
coating facility*

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Our quality criteria:

- German manufacturing expertise using only high-quality materials
- Hard, scratch-resistant precision anodizing
- Extremely narrow production tolerances

LINOS Nanobench

Like the Microbench, you can get this original from Qioptiq!

The Nanobench is a precise, extremely compact construction system for beam diameters of up to 1/2". Narrow production tolerances ensure that the optical axis is maintained and all components are mutually compatible.

With its large selection of optical elements, plus the option of integrating fiber couplers, the Nanobench has become one of the most popular construction systems in microoptics. You, too, can open up new applications with the Nanobench, and discover the highest precision with compact dimensions!

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Special Features:

Qioptiq offers a broad assortment of mounted optics for the Nanobench. Just look for the  symbol in the tables of optics.

Ideal Areas:

Fiber optics, microoptics, compact construction systems for applications with restrictive space requirements.

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Content

<i>Technical Features</i>	82
Nanobench Sets	
<i>Nanobench Mechanics Test Kit</i>	84
<i>Nanobench Set Mechanics</i>	85
<i>Nanobench Set Optics</i>	87
<i>Nanobench Basic Set</i>	88
Mounting Plates	
<i>Mounting Plate N</i>	89
<i>Mounting Plate N, latchable and rotatable</i>	90
Holder and Base Plates	
<i>Holder N 20</i>	90
<i>Holder N 40</i>	91
<i>Base Plate N</i>	91
<i>Base Plate N 40</i>	92
<i>Base Plate N 50</i>	92
Rods and Cubes	
<i>Set of Rods N</i>	93
<i>Cube N</i>	93
Positioners	
<i>X-Y Translation Stage N</i>	94
<i>Adjustment Plate N</i>	94
<i>Z-Adjuster N</i>	95
Fine Adjustment Units and Object Slides	
<i>Mirror Holder 45°</i>	95
<i>Mirror Holder 45°/7x12</i>	96
<i>Prism Holder</i>	96
<i>Rotary Support N</i>	97
Adapter	
<i>Microbench Adapter</i>	98
<i>Fiber Optic Adapter N FC</i>	98
<i>Fiber Optic Adapter N FSMA</i>	99
Tools and Screws	
<i>Tool Set N</i>	99
<i>Set of Set Screws M2.5x3</i>	100
<i>Set of Screws for Cube N</i>	100
<i>Set of Screws M2.5x6</i>	100
<i>Collar N</i>	100

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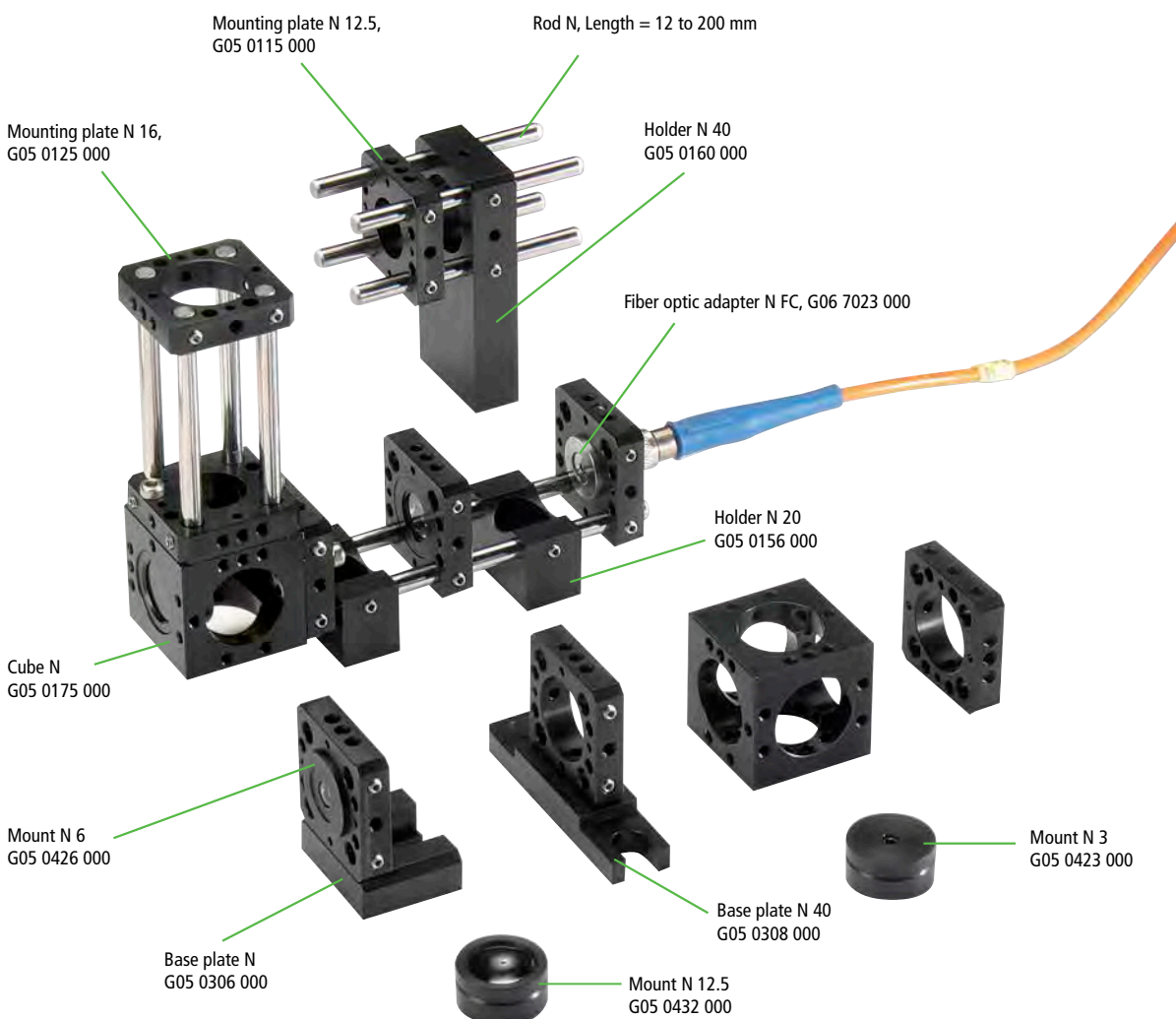
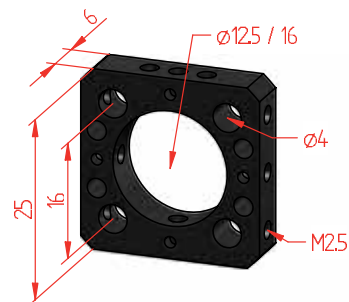
Technical Features

The Nanobench is a compact construction system for optics with diameters of up to 12.7 mm, and features exceptionally high precision. As with the Microbench system, mounted optics are fastened in mounting plates, which are connected by four rods to ensure highly precise alignment to the optical axis.

A variety of base plates makes it easy mounting Nanobench components directly on optical tables e.g. with its compact dimensions and flexibility in three-dimensional structures, the Nanobench is ideal for applications in micro-optics and fiber optics. Here, too, the LINOS catalog offers a broad selection of mounted optics that can be directly attached to the Nanobench. The  symbol helps to make it easy to find Nanobench-compatible products and accessories when ordering.

Adapters are available for fitting Microbench and Tube System C elements to our Nanobench system.

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- Ideal for use in micro- and fiber optics
- Compatible with the Microbench and Tube Mounting System C systems
- The ideal system for optics with \varnothing 3 mm to 12.7 mm
- Compact dimensions
- Modular design ensures greatest flexibility
- Extensive range of optics in Nanobench mounts

Basic construction elements

The functional principle of the Nanobench is identical to that of the Microbench system. The dimensions and tolerances, however, are on a smaller scale, compared to the Microbench's size. A wide range of mounting plates, holders, rods and cubes are available for creating sturdy and extremely compact systems.

Mounting plates

The bore holes for the rods on all Nanobench mounting plates have identical diameters and spacing. The edge length is 25 mm, and thickness is 6 mm. Where not indicated otherwise, the same dimensions apply for other components in the series.

- Lead-free aluminum alloy
- Matt black anodized surfaces
- Inner diameter: 12.5 or 16 mm with H7 tolerances

Four bore holes for mounting the connecting rods. The tolerances for the rod bore holes on the plates, together with the tolerances for the rods, guarantee a system with virtually zero play.

Rods

The rods are induction hardened to prevent marks made by adjustment screws. To ensure there is no play in the guidance of mounting plates and other components, the rods are lapped to fit.

- Hardened stainless steel with polished surfaces
- Lapped to fit
- Outer diameter: 4 mm
- Lengths: 12 mm to 200 mm

Cubes

Cube N is a multifunctional element for creating a beam diverter or integrating various optics in optic holders. Combined with rods and mounting plates, it lets you set up extremely compact three-dimensional structures. The mount diameter is 16 mm.

- Multifunctional elements for 3D structures
- Ideal for beam deflectors
- Inner diameter of bore holes: 16 mm
- M2.5 thread for mounting plates

The screws and threaded pins in the Nanobench system have M2.5 threads.

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Integration von Planoptik
Integration of planar optics

Mirror holder 45°/7x12
G05 0557 000
(with bonded beam-splitter plate 7x12 mm)

Mirror holder N G05 0568 000
(with bonded mirror 10x14 mm)

Rotation von Optik
Rotation of optics

Rotary support N in
Mounting plate N 16
with optics in mounts

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- For years now our customers are convinced by the quality of our bench systems. With this starter set we want to give you the chance to try the nanobench in your application.
- The Test Kit contains the basic components of the nanobench system such as rods and holders as well as all necessary screws and tools.
- In our webshop you will find all specific components for your application.

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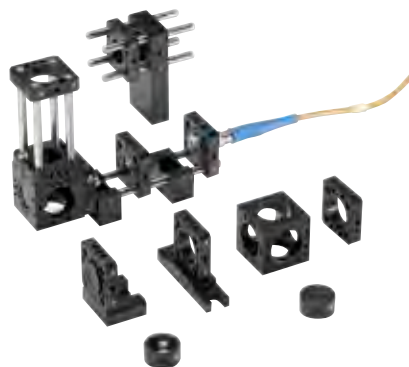
Please find further information about application examples, guidelines, frequently asked questions and more in our Brochure "Microbench, Bench and Rail systems - Introduction, solutions and applications". You will find the brochure for download at each Microbench product under "Docs + Drawings".

Mounting examples, the shown components are not identical to the content of the set

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- Includes all most important mechanic parts for an easy start with the Nanobench system.

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Nanobench Set Optics

Part No.	Qty.	Description
G05 2101 000	1	Plano-convex lens 2.5/3, mounted
G05 2102 000	1	Plano-convex lens 3/3, mounted
G05 2103 000	1	Plano-convex lens 5/5, mounted
G05 2105 000	1	Plano-convex lens 10/6, mounted
G05 2108 000	1	Plano-convex lens 15/10, mounted
G05 2109 000	1	Plano-convex lens 20/12,5, mounted
G05 2112 000	1	Plano-convex lens 50/12,7, mounted
G05 2114 000	1	Plano-convex lens 100/12,7, mounted
G05 2116 000	1	Plano-convex lens 200/12,7, mounted
G05 2201 000	1	Symmetric-convex lens 5/6, mounted
G05 2203 000	1	Symmetric-convex lens 10/10, mounted
G05 2206 000	1	Symmetric-convex lens 20/12,7, mounted
G05 2208 000	1	Symmetric-convex lens 50/12,7, mounted
G05 2210 000	1	Symmetric-convex lens 100/12,7, mounted
G05 2216 000	1	Best form lens, 6/3, mounted
G05 2217 000	1	Best form lens, 10/6, mounted
G05 2218 000	1	Best form lens, 20/10, mounted
G05 2302 000	1	Plano-concave lens, -6/6, mounted
G05 2303 000	1	Plano-concave lens, -10/6, mounted
G05 2305 000	1	Plano-concave lens, -10/10, mounted
G05 2308 000	1	Plano-concave lens, -50/12,7, mounted
G05 2309 000	1	Plano-concave lens, -100/12,7, mounted
G05 2402 000	1	Symmetric-concave lens, -20/12,7, mounted
G05 2403 000	1	Symmetric-concave lens, -50/12,7, mounted
G05 2404 000	1	Symmetric-concave lens, -100/12,7, mounted
G05 2003 000	1	Achromat, 6/3, mounted
G05 2004 000	1	Achromat, 10/6, mounted
G05 2006 000	1	Achromat, 20/10, mounted
G05 2007 000	1	Achromat, 25/12.5, mounted
G05 2010 000	1	Achromat, 40/12.5, mounted
G05 2012 000	1	Achromat, 50/12.5, mounted
G05 2014 000	1	Achromat, 80/12.5, mounted
G05 2015 000	1	Achromat, 100/12.5, mounted
G05 2018 000	1	Achromat, -10/6, mounted
G05 2017 000	1	Achromat, -20/8, mounted
G05 2019 000	1	Achromat, -35/12.5, mounted
G33 9911 000	2	90°-Prism, 10x10
G33 9952 000	1	90°-Prism, RAGV, 10x10
G33 5510 000	2	Beamsplitter cube, 10x10
G34 4140 000	2	Beamsplitter plate, 7x12
G34 0610 000	2	Plane mirror, ø10

- Includes all most important optic parts for an easy start with nanobench system
- Most optics already mounted for immediate use

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Nanobench Basic Set


It includes all components of the following cases:

*Nanobench Set Mechanics G050003000
Nanobench Set Optics G050004000*

- *Ideal set to start*
- *Most optics already mounted for immediate use*
- *Inclusive: Software-Package WinLens™ 3D (see chapter Optics Software for further information or on www.winlens.de)*

Nanobank-Grundkasten Nanobench Basic Set

Produktbezeichnung	Part No.
Item Title	Part No.
Nanobench Basic Set	G050006000

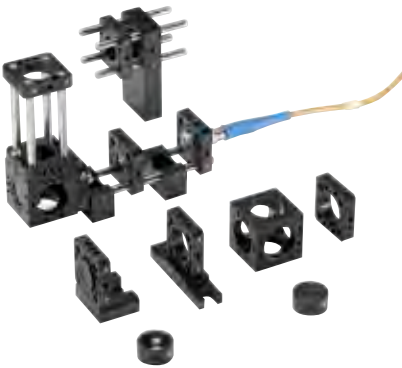
Mounting examples, the shown components are not identical to the contents of the set

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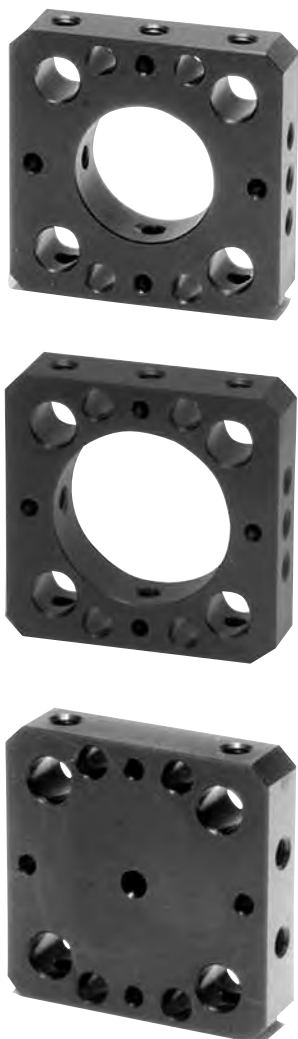
The Nanobench Basic Set is especially suited for a basic configuration of optical laboratories. It includes the most important optics and opto-mechanics as well as appropriate tools and screws. Additionally, a full version of our optical design software package WinLens™ is included. See more details in the chapter optic software or on www.winlens.de.

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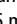
Mounting Plate N

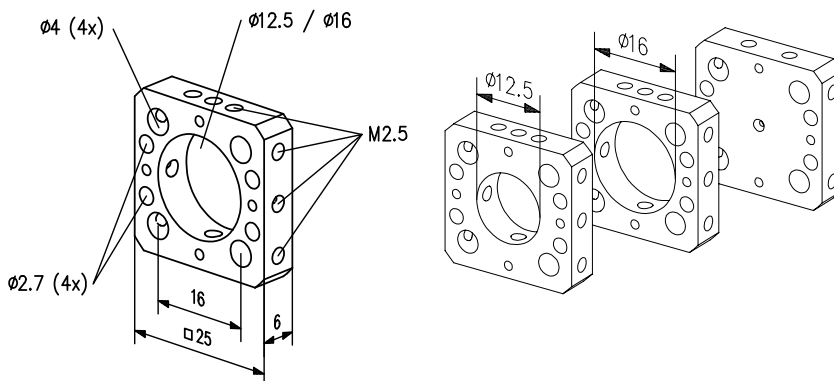- For 2, 3 and 4 rod setups
- To integrate mounts, adapters and connectors
- Available with tapped holes for attachment to Cube N
- For attaching rods and mounted lenses, you additionally require a set of Set screws M2.5x3 (G05 1520 000 or G05 1521 000), see tools and screws at the end of this chapter
- Set screws are included as standard components in Mounting Plate sets
- Mounted lenses for Mounting plate N 16 can be found in "Singlets" or "Achromats". All Nanobench mounts with an outer diameter of 16 mm are identified by an .

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

Produktbezeichnung	Part No.
Item Title	Part No.
Mounting Plate N 12.5	G050115000
Set of Mounting Plates N 12.5, 4 pcs.	G051115000
Mounting Plate N 16	G050125000
Set of Mounting Plates N 16, 4 pcs.	G051125000
Mounting Plate N without hole	G050111000

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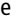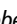
For compatible and appropriate optics (lenses, filters, etc) , see the Optics chapters. Optics compatible to our bench systems are marked as follows: Microbench  or Nanobench .

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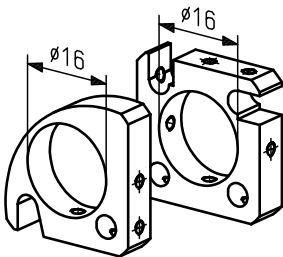Mounting Plate N, latchable and rotatable

- For later insertion into designs with rods
- Rotatable
- For attaching rods and mounted lenses, you additionally require a set of Set screws M2.5x3 (G05 1520 000 or G05 1521 000), see tools and screws at the end of this chapter

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
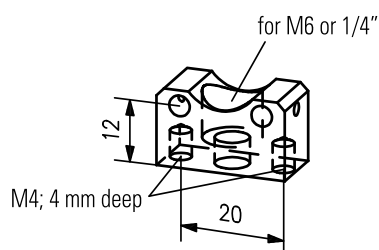
Holder N 20

- Holder for Nanobench setups
- To mount on plates or rods
- 20 mm optical axis height
- For attaching rods and mounted lenses, you additionally require a set of Set screws M2.5x3 (G05 1520 000 or G05 1521 000), see tools and screws at the end of this chapter
- Setscrews are included as standard components in Set of Holders N 20

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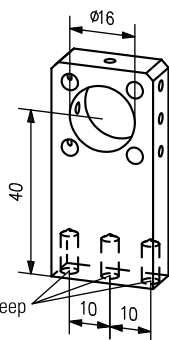
Holder N 40

- For insertion of mounts, adapters and connectors
- 40 mm optical axis height
- To mount on plates or rods
- For attaching rods and mounted lenses, a set of setscrews M2.5x3, G05 1520 000 or G05 1521 000 is required, see tools and screws at the end of this chapter
- Setscrews are included as standard components in Set of Holders N 40.
- Mounted lenses for Holder 40 can be found in "Singlets" or "Achromats". All Nanobench mounts with an outer diameter of 16 mm are identified by an .

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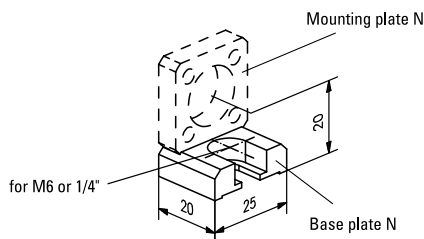
Base Plate N

- For attaching Mounting plate N to optical tables or posts
- 20 mm optical axis height
- Set of screws M2.5x6, G05 1523 000 is needed for fixing the Mounting plate N to the Base plate N, see Screws at the end of the chapter
- Screws are already included in Set of Base plates N.

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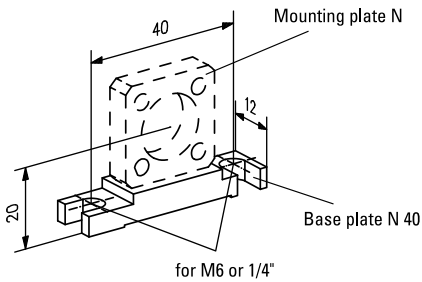
Base Plate N 40

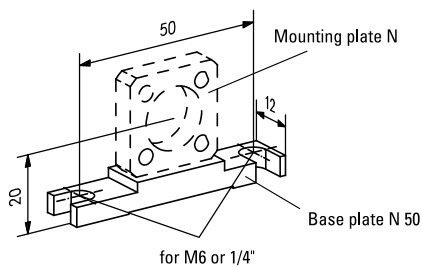- For fastening Mounting plates N onto Mounting plates with 40-mm tapped hole pattern
- 20-mm optical axis height
- Set of screws M2.5x6, G05 1523 000, for fixing the Mounting plate N to the Base plate N 40 required, see screws at the end of this chapter
- Screws are already included in Set of Base plates N 40.

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
Base Plate N 50

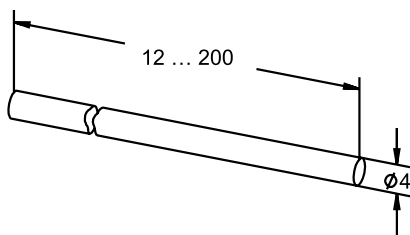- For attaching Mounting plates N onto Breadboards with 25-mm or 1"-tapped hole pattern
- 20-mm optical axis height
- Set of screws M2.5x6, G05 1523 000, for fixing the Mounting plate N to the Base plate N 50 required, see screws at the end of this chapter
- Screws are already included in Set of Base plates N 50.

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Set of Rods N

- For connecting different components such as Mounting Plate N or Cube N
- To realize 2, 3, and 4 rod setups
- Acid resistant stainless steel
- Induction hardened, lap-ground surfaces

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
Cube N

- For assembling compact and rigid 3-D setups
- Useful for creating right angles in optical setup
- For insertion of beamsplitters, mirrors, and prisms you will need Mirror Holder 45° or the Prism Holder.
- For attaching rods you will need Mounting Plate N.
- Set of screws (G05 1522 000) is included.

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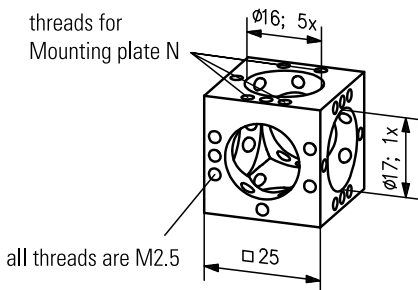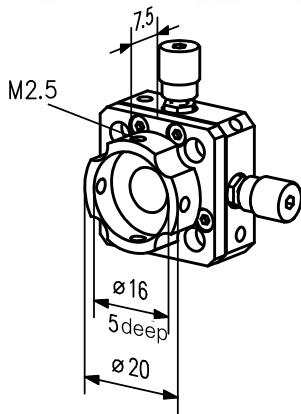
For compatible and appropriate optics (lenses, filters, etc) , see the Optics chapters. Optics compatible to our bench systems are marked as follows: Microbench or Nanobench .

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X-Y Translation Stage N

- Mounting diameter 16 mm
- 8 mm aperture
- Adjustment screws with hex socket head screws
- Fine adjustment screws with 0.25 mm pitch
- ± 0.5 mm alignment range

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The XY translation stage N is ideally suited for fiber coupling applications. Please find compatible adapters in the adapter section of this chapter.

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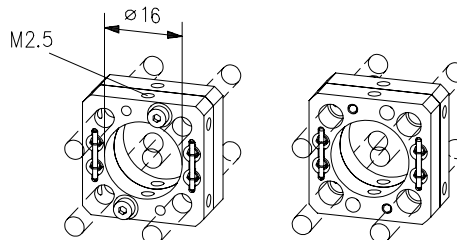
Adjustment Plate N

- Two plates connected with springs.
- Manual adjustment in X and Y direction
- Lockable

- Mounting diameter of 16 mm
- ± 0.5 mm alignment range

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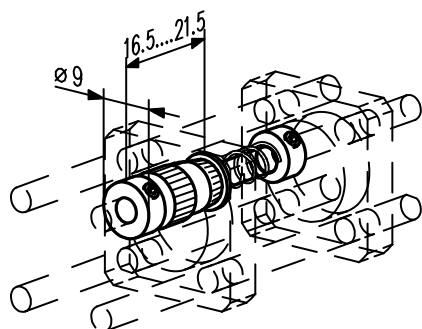
Z-Adjuster N

- Precision translation unit with 5 mm travel
- 0.25 mm pitch
- Backlash free adjustment of optics, pinholes, fiber connectors etc.
- Compact design
- Can be inserted into existing setups

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
- Consisting of a micrometer, compression spring and collar

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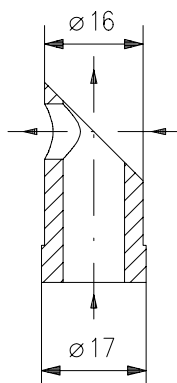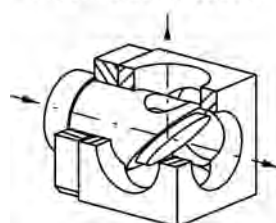
Mirror Holder 45°

- Accepts bonded elliptical mirrors or beamsplitter plates, e.g. 12.5x18 mm
- Insert in Cube N or Mounting plate N 16
- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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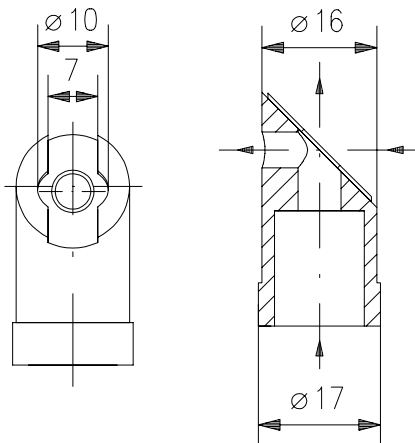
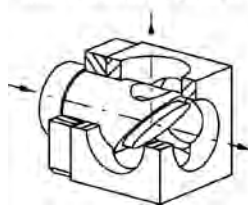
Mirror holder in Cube N


Mirror Holder 45°/7x12

- Accepts bonded rectangular beamsplitter plates 7x12 mm
- Countersunk for circular mirrors with 10 mm outer diameter
- Insert in Cube N or Mounting plate N 16
- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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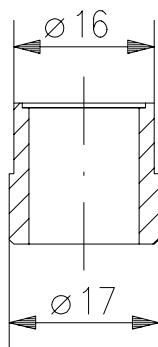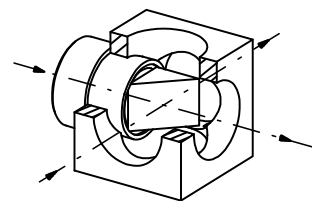
Prism holder in Cube N


Prism Holder

- To bond prisms, beamsplitter cubes or thin film polarizers with 10x10 mm dimensions
- To insert in Cube N or Mounting plate N 16
- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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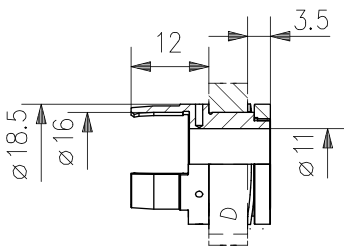Rotary Support N

The Rotary support is designed to be clipped into Mounting Plate N 16 and is tensioned by a disc spring to ensure backlash-free rotation. A lateral bore hole ($\varnothing 1.1$ mm) is provided for insertion of an optional lever to facilitate turning the support. Rotary Support N can also be used as a quick-change set for optics in 16 mm mounts, as the mounted optics are simply clipped into place.

- Insert for mounting plates N 16
- Easy plug in of optics in mount N 16
- Free from backlash by use of a spring
- Also usable for quick exchange of different optics in mount N 16
- Rotatable by lever (not included) in sideways bores with $\varnothing 1.1$ mm
- Mounting plate N16 and optics not included

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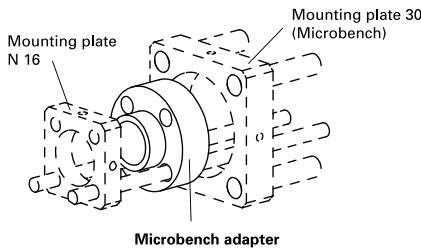
Microbench Adapter

- To connect the LINOS cage mounting systems Microbench and Nanobench with Mounting plate N 16 and Mounting plate 30 (Microbench)
- Through hole for Nanobench rods

Mikrobankadapter Microbench Adapter

Produktbezeichnung	Außen-Ø *) (mm)	Innendurchmesser (mm)	Länge (mm)	Part No.
Item Title	Outer-Ø * (mm)	Inner diameter (mm)	Length (mm)	Part No.
Microbench Adapter	25	16	10	G050305000

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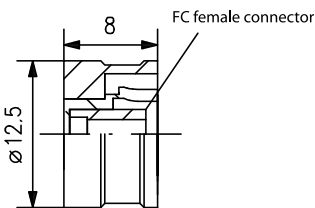
Fiber Optic Adapter N FC

- Mounting flange without optics
- To connect fibers with FC connector
- Outer diameter 12.5 mm to insert into Mounting plate N 12.5
- Outer diameter 16 mm to insert into Mounting plate N 16 or XY Adjustment plates N

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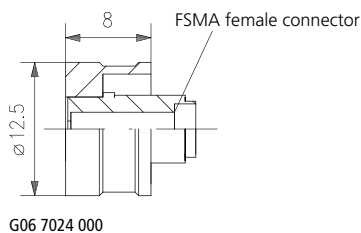
Fiber Optic Adapter N FSMA

- Mounting flange without optics
- To connect fibers with FSMA connector
- Outer diameter 12.5 mm to insert into Mounting plate N 12.5
- Outer diameter 16 mm to insert into Mounting plate N 16 or XY Adjustment plates N

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Tool Set N

- Consists of:
- Tube wrench for retaining ring M13.5
 - Allen wrench, size 1.3 mm
 - Allen wrench, size 2 mm
 - Tweezers
 - Allen wrench with grip and spherical head, size 1.3 mm
 - Allen wrench with grip and spherical head, size 2 mm

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Produktbezeichnung	Part No.
Item Title	Part No.
Tool Set N	G050501000

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Set of Screws for Cube N

For attaching Mounting plates N to Cubes N
Consists of:

- Setscrews M2.5x3, 10 ea.
- Hexagon socket head cap screws M2.5x8, 10 ea. (DIN 912)

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Set of Screws M2.5x6

- For fastening Mounting plates N to Base plates N
- Quantity 25

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Collar N

- For use as clamps or adjustable stops

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Our quality criteria:

- *German manufacturing expertise using only high-quality materials*
- *Hard, scratch-resistant precision anodizing*
- *Extremely narrow production tolerances*

LINOS Tube System C

In addition to the open Microbench system, Qioptiq also offers a closed, modular Tube Mounting System.

The LINOS Tube Mounting System C is a component construction system for fast and trouble-free set-up of laboratory instruments, prototypes and small batch series. Tubes are available in diameters of 30 mm and 35 mm. Threaded sockets and threaded rings enable connection of a wide range of optics, with diameters from 6 mm to 31.5 mm. With the Tube Mounting System C, you can create your own lens and expander systems.

Adapter rings are available for compatibility with the connecting threads on C-mounts, microscopes and CCD cameras. The C30 Tubes feature C-mount connection threads and are particularly well suited for setting up individual CCD camera lenses.

Fine adjustments and focusing mounts are available to enable minimal length adjustments for high-precision focusing. The LINOS Tube Mounting System and the LINOS Microbench are excellent in combination; for example, to create a light-tight and dust-tight beam path.

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Special Features:

Mount system for our optics with diameters of up to 31.5 mm, for easy assembly of existing lenses and fast setup of prototypes. Continuous system expansion by adding new components.

Ideal Areas:

Light-tight and dust-tight beam paths; rapid setup of existing lenses; prototypes; small batch series.

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Content

<i>Technical Features</i>	104
Set	
<i>Mechanics Set</i>	106
Tubes C 30	
<i>Tube C30</i>	107
<i>Closing Tube C30x6</i>	107
Tubes C 35	
<i>Tube C35</i>	108
<i>Lens Tube C35x50</i>	108
<i>Retaining Mount 31.5 mm</i>	109
Cubes and Cube Inserts	
<i>Beam Deflector Cube C</i>	109
<i>Beam Deflector Cube C-5</i>	110
<i>Lid C</i>	110
<i>Beamsteering Mirror Mount, adjustable</i>	111
<i>Prism Table Insert, Adjustable</i>	111
<i>Beamsplitter Cube Insert, Adjustable</i>	112
Mounts and Connecting Elements	
<i>Spatial Filter Module C30x37,5</i>	113
<i>Filter Support C</i>	114
<i>Threaded Mounts</i>	115
<i>Iris Diaphragm Set</i>	116
<i>Light Trap</i>	116
<i>Beam Dump</i>	116
<i>Optic Mount 1", C30x15</i>	117
<i>Mount 31.5</i>	117
<i>Adjusting Tube C30 with Locking Collar</i>	118
<i>Focusing Mount C30</i>	118
<i>Connecting Tube C30 FC / C30 FSMA</i>	119
<i>Fine Adjustment C30</i>	119
<i>Focusing Mount C35</i>	120
<i>Connecting Tube C30 with Adjustable Mount 6</i>	120
<i>Eyepiece Adapter C30</i>	120
<i>Tube C with DS Polarizer, Rotatable</i>	121
<i>C-Mount Adapter M17</i>	121
Retaining Rings, Spacers and Tools	
<i>Adapter 1" to W0.8" or Ø 25 mm</i>	122
<i>Adapter W0.8" to 1"</i>	122
<i>Knurled Ring C30x8</i>	122
<i>C-Mount Microbench Adapter</i>	123
<i>Adapter Ring C</i>	123
<i>Adapter Ring C-W0.8</i>	123
<i>Retaining Ring M32.5x0.75</i>	124
<i>Retaining Ring M32.5x0.75-8</i>	124
<i>Spacer Ø 31.5</i>	124
<i>Retaining Ring M23.2x0.75</i>	124
<i>Spacer Ø 22.4</i>	125
<i>Flat Wrench</i>	125
<i>Tube Wrench 22.4/31.5</i>	125
Clamp Holder	
<i>Tube C - Clamp Holder</i>	126

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Combination of Tube Mounting System C and Microbench

Optics adjustment along the Z axis

Adapter C-Mount/Microbench
G06 1659 000

Tubes C30 with optics
Focusing mount C35,
G06 5239 000

Technical Features

The Tube Mounting System C is a modular component construction system for fast and trouble-free set-up of laboratory instruments, prototypes and small-batch production. Tubes are available in diameters of 30 mm and 35 mm. Threaded sockets and threaded rings enable the connection of a wide range of optics, with diameters from 6 mm to 31.5 mm.

With the Tube Mounting System C, creating your own lens and expander systems is a breeze. Adapter rings make this system compatible with the connecting threads on C-Mounts, microscopes and CCD cameras. Fine adjusters and focusing mounts let you make small changes in the length for highly precise focusing. Tube Mounting System C and Microbench are excellent in combination; for example, to create a light-tight and dust-tight beam path.

The Tube Mounting System C is compatible to the Microbench and is being constantly extended with new components.

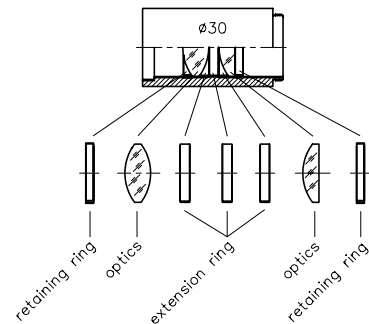
- To realize your laboratory instruments, samples, prototypes and small-batch production, including lenses and expander systems
- Modular component system
- For unmounted optics with \varnothing 6 mm to 31.5 mm
- Compatible with C-Mount, microscope and CCD camera threads

Examples of Optics Assemblies

Lens System in Tube Mounting System \varnothing 35 mm


Lens System in Tube Mounting System \varnothing 30 mm


Examples of Complete Systems

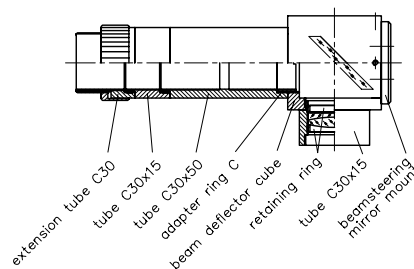
CCD-Camera Objective with Bent Beam Path

Beam Expander System

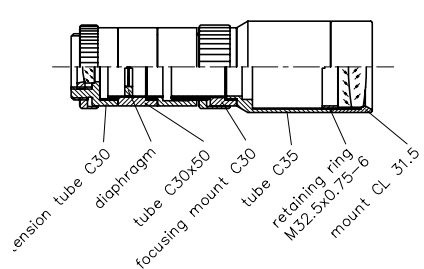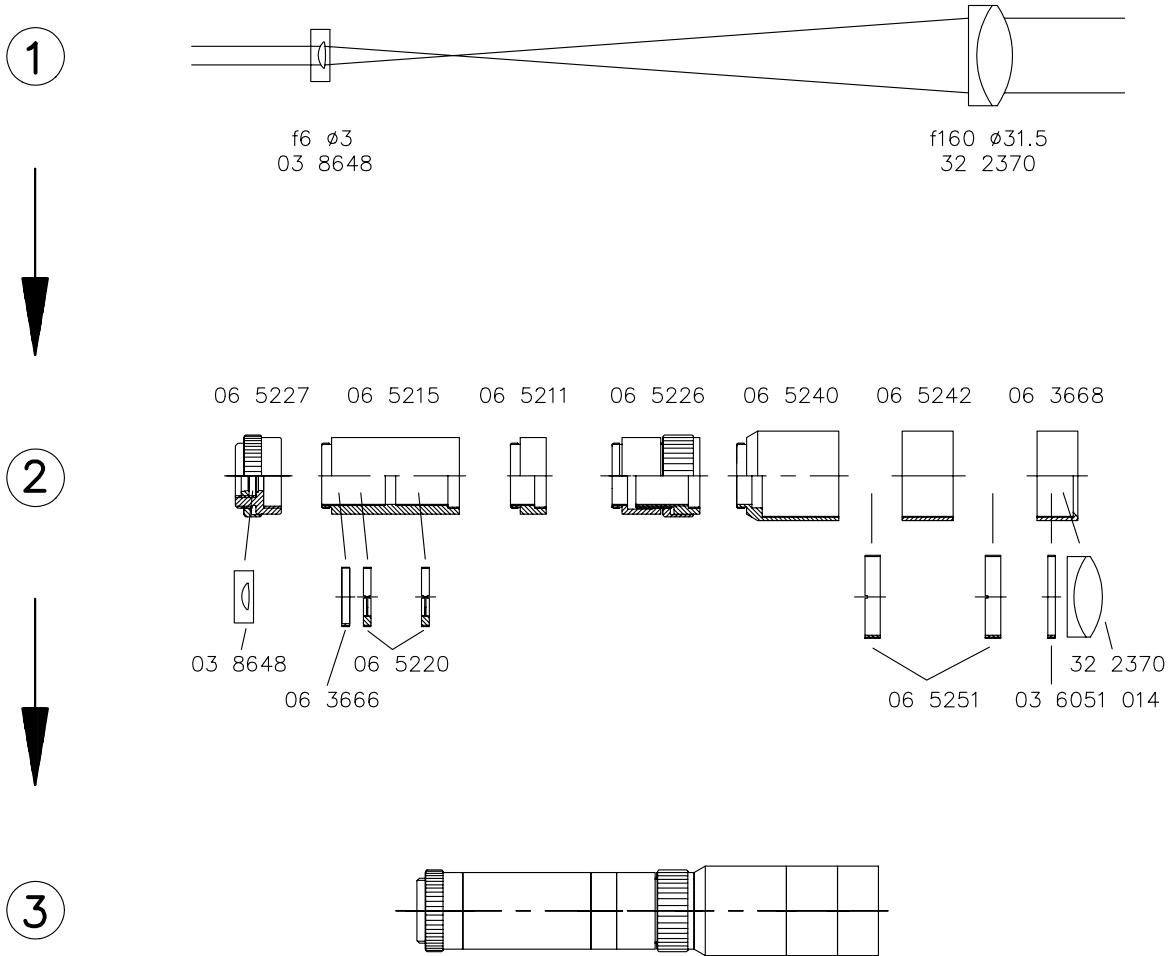Example: Construction of a Laser beam expander system 26x with components of the Tube Mounting System C

Specifications:

Expansion ratio	26x		
Entrance optic	f=6 mm	Ø 3 mm	free Ø 2.5 mm
Exit optic	f=160 mm	Ø 31.5 mm	free Ø 30 mm
Wavelength	450-700 nm		
Length	180 mm		
Diameter	35 mm		
focusable, with locking collar			
Mounting thread	1-32UN (C-Mount)		

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Tube Mounting System C - Mechanics Set

part No.	Qty.	Description
G065211000	2	Tube C30x10
G065212000	3	Tube C30x15
G065215000	1	Tube C30x50
G065218000	1	Tube C30x85
G061652000	1	Iris diaphragm C15/30
G065226000	1	Focusing mount C30
G065225000	1	Adjusting tube C30 with locking collar
G065230000	1	Eyepiece adapter C30
G065220000	1	Iris diaphragm set
G065235000	1	Beam deflector cube C
G061667000	1	Adapter ring C
G065096000	1	Beamsteering mirror mount, adjustable
G065097000	1	Prism table insert, adjustable
G065098000	1	Beamsplitter cube insert, adjustable
G065240000	1	Tube C35
G065242000	2	Tube C35x20
G065244000	1	Tube C35x30
G061653000	1	Iris diaphragm C22/35
G063668000	2	Mount CL 31.5
G036051014	5	Retaining ring M32.5x0.75
G065251000	5	Retaining ring M32.5x0.75-6
G066051013	5	Extension ring Ø31.5
G063666000	10	Retaining ring M23.2x0.75
G063010012	10	Extension ring Ø22.4
G061097000	1	Flat wrench 22.4
G061099000	1	Flat wrench 31.5
G061090000	1	Screw driver 1.8

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- Includes all most important mechanic parts for an easy start with the Tube mounting system.

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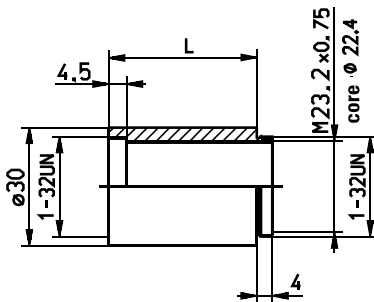
Tube C30

- To mount optics with outer Ø 22.4 mm with two retaining rings M23.2x0.75 (G06 3666 000)
- With threaded mounts lenses and achromats 6 to 19 mm Ø can be mounted into the tubes
- External-Ø 30 mm
- Internal and external 1-32UN (C-Mount) thread
- Internal thread M23.2 x 0.75 for optics mounting (for L = 85 mm: on both sides internal thread cut of 30 mm depth)
- Internal-Ø 22.4 mm

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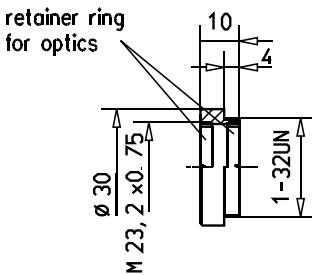
The adaption thread of the tube system C is the so-called "C-mount", a standard thread for camera objectives. The diameter is 1", the pitch 1/32". The standard distance between image plane and flange of the objective thread is 0.69" (17.526 mm).

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Closing Tube C30x6

- Mount for filters, protective disks, auxiliary close-up lenses etc.
- End ring for Tube C components (to cover C-Mount)
- Two retaining rings for optics with 22.4 mm diameter are included

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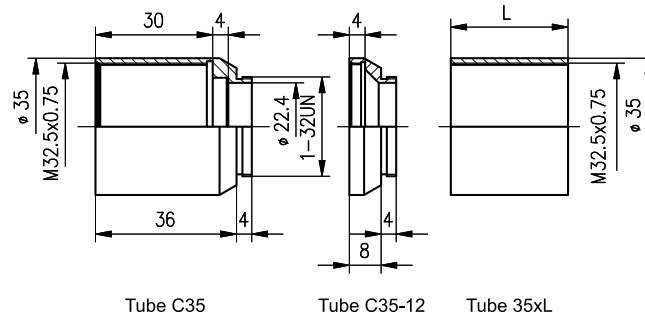
Tube C35


- Outer-Ø 35 mm
- Tube C30 with 1-32UN (C-Mount) thread
- Internal thread M32.5x0.75
- Distance fixing via spacer (available with 2mm, 5mm and 10mm width)
- Use suitable retaining- and extension rings for Ø 31.5 mm pinholes, reticles and optics
- Retaining ring (G06 5251 000) is necessary for coupling

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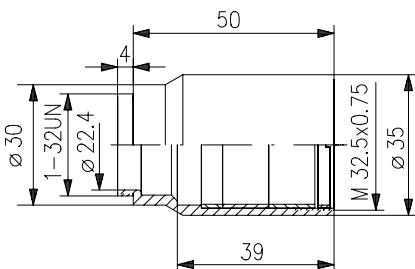Lens Tube C35x50


- For assembling simple lenses
- With retaining ring M32.5x0.75
- Extension rings with 5mm and 10mm width are included

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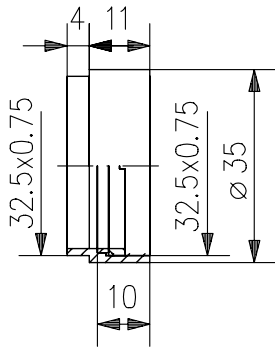Retaining Mount 31.5 mm

- For filters, polarization filters, dispersion plates, etc. with Ø 31.5mm
- For all tube elements with M32.5x0.75 thread
- Stackable thanks to mounting thread on both ends
- Complete with retaining ring and premounting ring

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
Beam Deflector Cube C

- For beamsteering mirror mount and prism table insert
- 40x40x40 mm, rounded corners
- 2 x internal thread C (1-32UN)
- 1 x Ø 30 mm
- Equipped with 1 adapter ring C (G06 1667 000)
- Use Connecting tube (G06 5225 000) or Focusing mount (G06 5226 000) for optimum alignment when combining two cubes.

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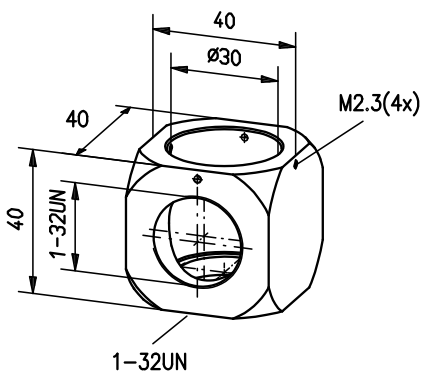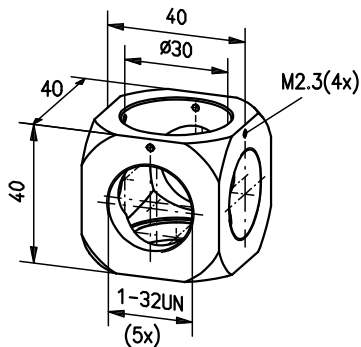
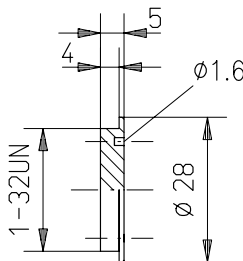
Beam Deflector Cube C-5

- For beamsteering mirror mount and prism table insert
- 40 x 40 x 40 mm, rounded corners
- 1 x Ø 30 mm
- 5 apertures with internal thread C (1-32UN)
- Equipped with 1 adapter ring C, G06 1667 000
- 3 lids, G06 5200 000, included

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
Lid C

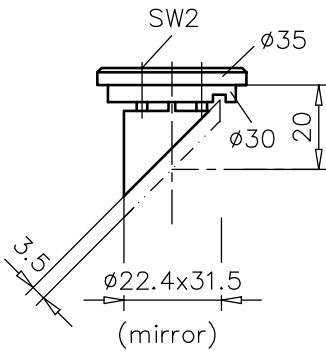- To cover CCD cameras, tubes, cube apertures etc. with internal thread C (1-32UN)

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

Beamsteering Mirror Mount, adjustable

- 90° beam deflection
- For mirror with Ø 22.4 x 31.5 mm e.g. G340003000
- Can be adjusted by spring joint
- Can be used in cube C
- Compatible with Microbench system via Ø 30 mm

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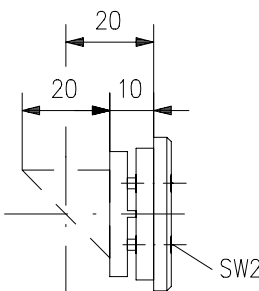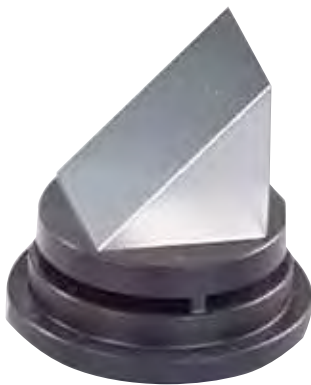Prism Table Insert, Adjustable

- To insert 90° prisms and splitter cubes with 20 mm edge length
- 1 planar side for gluing mirrors or plates
- Adjustable via spring joint
- You will find prisms, e.g. G339953000, in chapter plano optics, section prisms
- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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
Beamsplitter Cube Insert, Adjustable

- For beamsplitter cubes 20 x 20 x 20 mm
- Adjustable with spring joint
- To insert in cube C
- You will find beamsplitter cubes, e.g. G33 5520 000, in chapter plano optics, section beamsplitters.
- For instructions on gluing your optics onto the holder, please see section "Docs + Drawings". We are happy to take on this task for you. Please contact your local sales partner and we will provide an individual quote.

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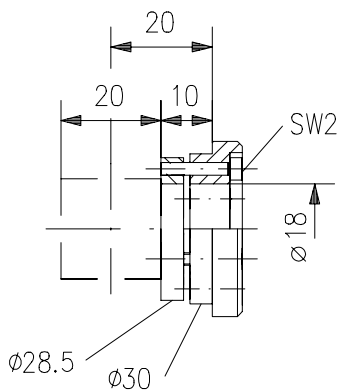Please find further information about application examples, guidelines, frequently asked questions and more in our brochure "Microbench, Bench and Rail systems - Introduction, solutions and applications". You will find the brochure for download at each Microbench product under "Docs + Drawings".

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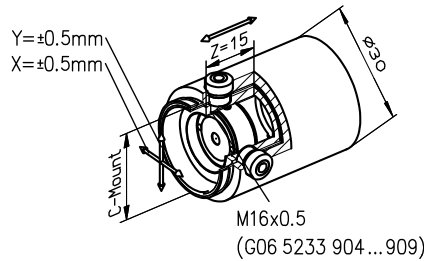
Spatial Filter Module Tube C30x37.5

The spatial filter module provides the possibility to set up beamexpanders, with spatial filters, with the tube mounting system. It provides precise xy-adjustment as well as adjustment along the optical axis (z).

- For spatial filter use
- C-mount thread on both sides
- Inner thread M16 for mounted pinholes
- ± 0.5 mm travel
- Spring loaded, backlash free adjusters

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Filter support including Filter holder for 1 filter

Filter support including Filter holder for 3 filters

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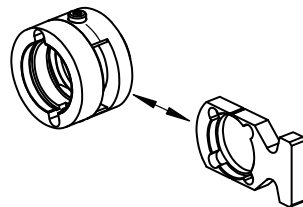
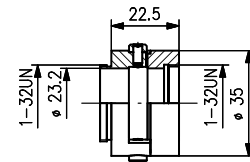
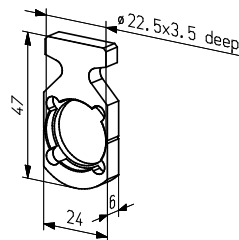
Filter support, including Filter holder for 1 filter, alternatively Filter holder for 3 filters


Filter holder for 1 filter, accessory to the Filter support

Filter Support C

Filter support for fast exchange between different filters, without disassembling the tube system. The filter will be glued into the holder and can be slid into the support.

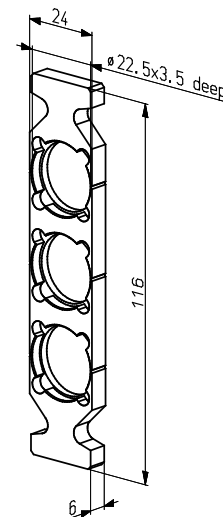- For unmouted filter diameter 22.4 mm.
- Holder in version for 1 or 3 filters available

Filter holder for 3 filters, accessory to the Filter support

Filterschieber C

Filter Support C

Produktbezeichnung	Part No.
Item Title	Part No.
Filter Support including Filter holder for 1 filter	G065255901
Filter Support including Filter holder for 3 filters	G065256901
Accessory, Filter Holder for 1 filter	G065255000
Accessory, Filter Holder for 3 filters	G065256000

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
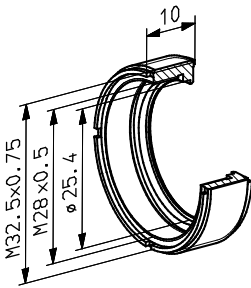
Threaded Mounts

- To screw in Tube C30 respectively Tube C35
- External M23.2 x 0.75 thread or M32.5 x 0.75 (G063 698 000)
- For optics with diameter 6 to 19 mm or diameter 25.4 mm (G06 3698 000)
- Diameter 19 mm (G063 689 000) for reticles
- Threaded Mount G063 689 000 also suitable for reticles with Ø 19 mm
- See the "Singlets" and "Achromats and Lens Systems" sections for details on a variety of unmounted optics.
- For reticles, see "Reticles/Test Targets" under "Plano Optics."

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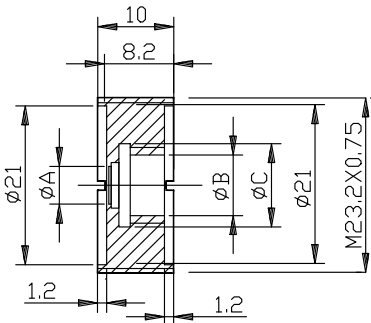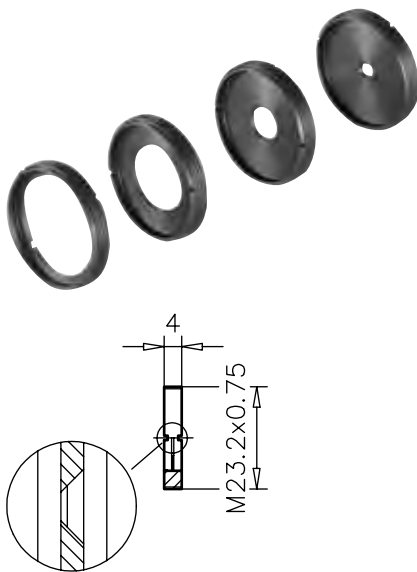
The tube system C is ideally suited to build individual beam expanding systems. An example is given in our brochure "Microbench, Bench and Rail systems". You will find the brochure for download at each Microbench product under "Docs + Drawings".

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
Iris Diaphragm Set

Set of 6 diaphragms with \varnothing 2.5 - 4 - 6 - 9 - 12 - 18 mm

- For use as stops or light traps
- M23.2x0.75 thread
- Suitable wrenchs Tube wrench (G06 1100 000) or Flat wrench 22.4 / 31.5 (G0 61097 000) not included.

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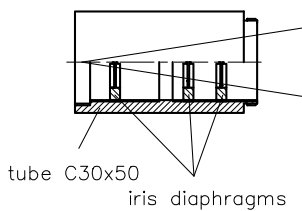Beam Dump

The beam dump is a solid aluminum cone, with a fluted, black anodized surface. The length is 46 mm. The cone can be mounted into components of the Tube Mounting System with the thread M23.2x0.75.

In conjunction with a tube C30x50 (G065215000) a beam dump for low power can be built.

- Fluted, black anodized surface
- Massive aluminum
- Connection thread M23.2x0.75

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

Optic Mount 1", C30x15

- Mount for 1"- optic
- Equipped with retaining and extension ring
- Two lenses can be mounted one by one (extension ring width 5 mm)

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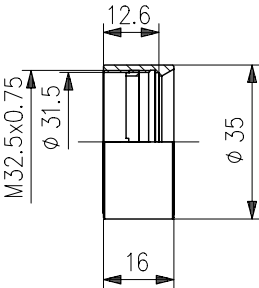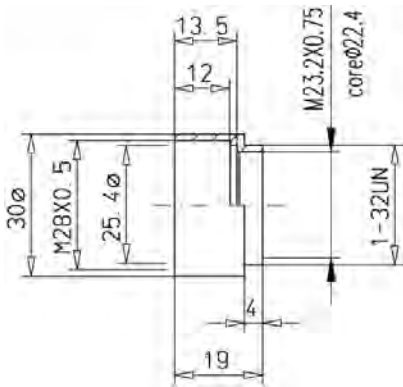Mount 31.5

- For Ø 31.5 mm optics
- Equipped with retaining
- For stacking two Mounts extension ring G065251000 is required

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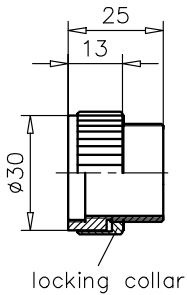
CAD files of all mechanical components can be found on the respective product sheets at www.qioptiq.com below the button "Docs and Drawings" .

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
Adjusting Tube C30 with Locking Collar

- Suitable for CCD cameras
- For connecting two cubes
- Additional mounting of optics with $\varnothing 22.4$ possible
- Easy locking of position by locking collar

- Tube- \varnothing 30 mm
- Tube length 25 mm
- External 1-32UN (C-Mount) thread, 15 mm long with locking collar \varnothing 32 mm
- Travel max. 10 mm
- Internal 1-32UN (C-Mount) thread
- Inner- \varnothing 22.4 mm with thread M23.2x0.75 for mounting additional optics

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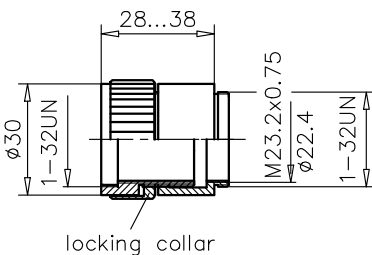Focusing Mount C30

- Like adjusting tube C30, but with additional Tube C30x15
- Inner- \varnothing 22.4 mm with thread M23.2x0.75 for mounting additional optics
- For connecting two cubes
- Additional mounting of optics with $\varnothing 22.4$ possible
- Easy locking of position by locking collar

- Tube- \varnothing 30 mm
- Adjustable tube length 28 - 38 mm
- External connection 1-32UN (C-Mount) thread
- With locking collar, outer- \varnothing 32 mm
- Internal connection 1-32UN (C-Mount) thread
- Inner- \varnothing 22.4 mm with thread M23.2x0.75 for mounting additional optics

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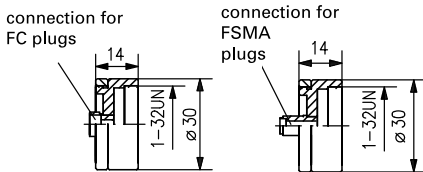
Connecting Tube C30 FC / C30 FSMA

Mounting flange without lens

- For connection of fiber light guides with FC/FSMA plugs
- C-Mount threads on both ends (1-32UN)
- Male threads covered by sleeve nuts
- Outer Ø 30 mm

Anschlussstube C30 FC / C30 FSMA Connecting Tube C30 FC / C30 FSMA

Produktbezeichnung	Part No.
Item Title	Part No.
Connecting Tube C30 FC	G067056000
Connecting Tube C30 FSMA	G067057000

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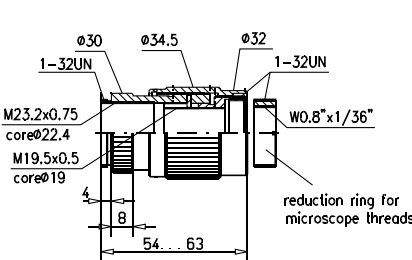
Fine Adjustment C30

- For z-axis adjustment of optics
- For connection of microscope lenses with adapter G06 1668 000 (included)
- Travel 3 mm or 6 mm
- Internal 1-32UN thread
- External 1-32UN thread

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
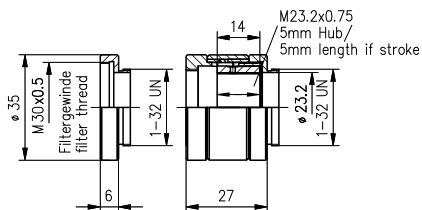
Focusing Mount C35

- For linear translation of optics in threaded mounts with thread M23.2x0.75
 - Direct mounting of optics with 22.4 mm diameter between two retaining rings
 - Length of the part stays same while focusing
 - Filter adapter included
 - Translation can be fixed with setscrew
- C-Mount thread on both sides

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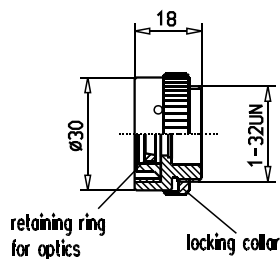Connecting Tube C30 with Adjustable Mount 6

- Adjustable mount of optics with 6 mm diameter
 - Application: For entrance optics of expanding systems or fiber light guides
- Mounting with included retaining ring
 - Centering range ± 0.5 mm

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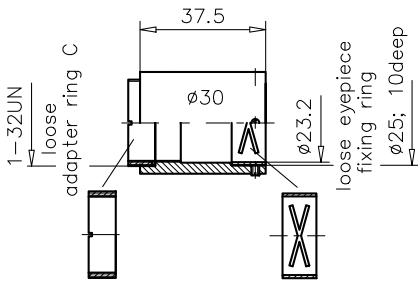
Eyepiece Adapter C30

- Outer-Ø 30 mm
- Inner-Ø 22.4 mm / Ø 25 mm with insert for Ø 23.2 mm eyepieces
- Loose adapter ring with thread C for either internal or external thread

Okularadapter C30
Eyepiece Adapter C30

Produktbezeichnung	Part No.
Item Title	Part No.
Eyeiece Adapter C30	G065230000

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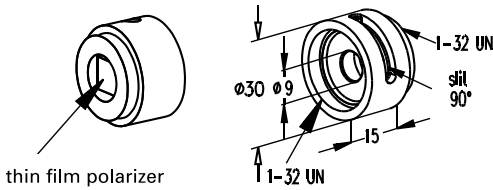
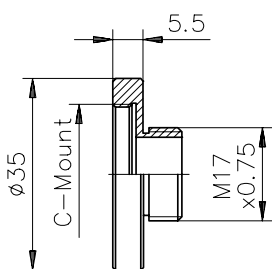
C-Mount Adapter M17

- For fitting objectives to C-Mount threads
- For connecting objective lenses with mounting thread M17x0.75

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
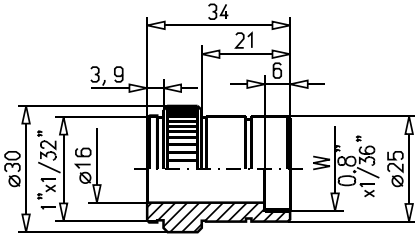
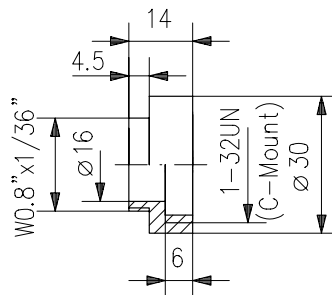
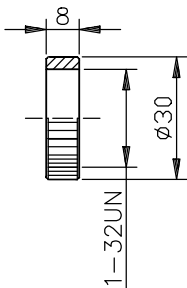
Knurled Ring C30x8

- For connecting two male threads
- Internal thread C (1-32UN)
- To change from male to female thread

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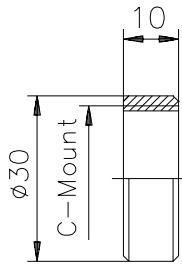
C-Mount Microbench Adapter

- Fits components with C-Mount thread to Ø 30 Microbench mounting plates

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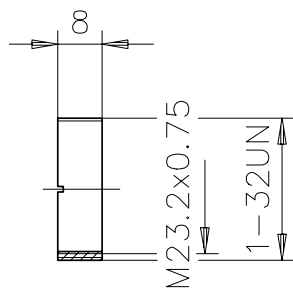
Adapter Ring C

- For connecting 2 threads C
- External thread C (1-32UN) 8 mm wide
- Internal thread M23.2x0.75 with inside-Ø 22.4 mm

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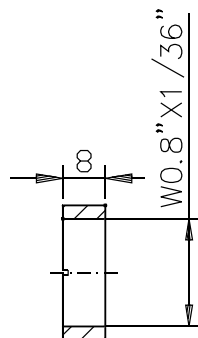
Adapter Ring C-W0.8

- To connect microscope objectives to tubes
- External thread C-Mount (1-32UN)
- Internal thread W0.8"
- 8 mm wide

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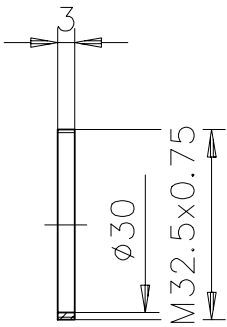
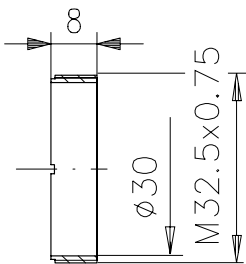
Retaining Ring M32.5x0.75-8

- To connect two $\varnothing 35$ mm tubes
- To connect $\varnothing 35$ mm tube to Mount CL 31.5 (G06 3668 000)

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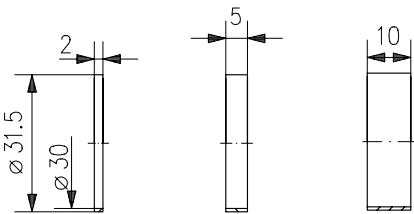
Spacer $\varnothing 31.5$

- For distance fixing of $\varnothing 31.5$ mm optics
- Different types with 2, 5 and 10 mm width

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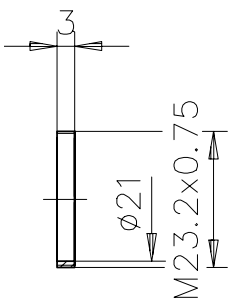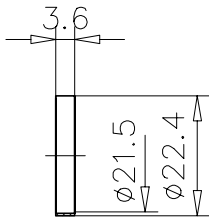
Flat Wrench

- For mounting retaining rings and diaphragms
- For Ø 22.4 mm or 31.5 mm

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Tube Wrench 22.4/31.5

- For mounting components with Ø 22.4 / 31.5 mm
- Suitable for retaining rings M23.2 x 0.75 and M32.5 x 0.75

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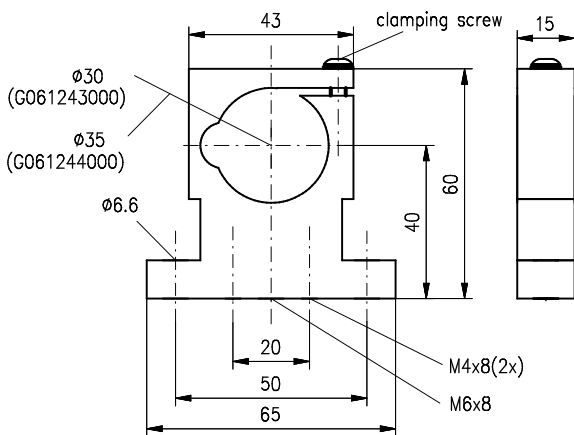
Tube C - Clamp Holder

- To mount components with outer \varnothing 30 or 35 mm
- Universal bores for mounting on Microbench mounting plates, breadboards or optical benches
- 40 mm centerline height
- Inner \varnothing 30 or 35 mm

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Profit from our 130 years of experience in the production of precision optical and mechanical components.

Früher / In former times

Montage und Mechanikfertigung in Göttingen um 1910
Mounting and production of mechanical components in Goettingen about 1910

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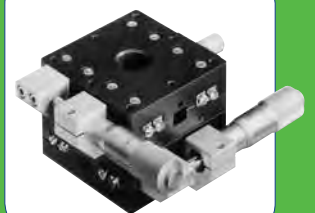
Our quality criteria:

- Custom production
- Precise micrometers
- Use of high-quality guides with maximum evenness

LINOS Positioners

Are you trying to find a solution for your positioning requirements? You can find it at Qioptiq!

We have a broad range of positioner types, including manual linear positioners, X-Y stages and rotary stages with assorted adjustment ranges and load capacities.

The manual LINOS positioners are equipped with micrometer screws with 0.25 or 0.5 thread pitches, which permit an adjustment sensitivity of better than 1 µm.

Piezosystems with suitable controllers complement the portfolio of our positioning systems.

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Special Features:

In addition to a large selection of sizes, adjustment ranges and load capacities, LINOS also offers the option of replacing the micrometers with actuators.

Ideal Areas:

Precision positioning of optics; microscopy.

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Content

<i>Technical information</i>	132
Manual Positioners	
<i>Linear Positioners TM / TMS</i>	134
<i>Linear Stages TB</i>	136
<i>Mounting Bracket WM 50-1/80-1 for Linear Stages TB</i>	138
<i>Mounting Bracket WM 50-2/80-2 for Linear Stages TB</i>	139
<i>Linear Stages T 10 M/F and TS 10 M/F</i>	140
<i>Linear Stages TL 45 / TL 65</i>	142
<i>Linear Stages TL 45 / TL 65 with Digital Micrometer</i>	143
<i>Linear Stages TL 65 with Side Drive</i>	144
<i>Mounting Components</i>	145
<i>X-Y Stages XY 60 / XY 85</i>	146
<i>X-Y Stage XY 85-25 with Digital Micrometer</i>	147
<i>X-Y Stages XY 85 with Side Drive</i>	148
<i>X-Y Stages XY 200</i>	149
<i>X-Y Stages XY 200 with Digital Micrometer</i>	150
<i>Clamping Set for Stages</i>	151
<i>Goniometer Stages 60-60 / 60-80</i>	152
<i>Rotary Stages RT 60 / RT 90 / RT 120</i>	153
<i>Z-Axis Stages HT 60-5 / 90-5 / 120-10</i>	154
<i>Vertical Translation Stage HT 100-40</i>	155
<i>Tilting Stage NT 60-3</i>	156
<i>Prism Stages 40 S / 60 S</i>	157
Micrometers and Adjustment Screws	
<i>Micrometers G</i>	158
<i>Precision Adjustment Screws G, bushing with connection thread</i>	158
<i>Fine-Adjustment Screws, Flat Bushing</i>	159
<i>Precision Lead Screws Lees</i>	160
Piezo Systems	
<i>CN 30 Controller for Piezo Electric Inertial Motors</i>	161
<i>CF 30 Controller for Piezo Electric Inertial Motors</i>	161
<i>Connection cables for piezo positioners</i>	161
<i>MS 30 Miniature Translation Stage</i>	162
<i>MS 38 Miniature Translation Stages with Free Opening</i>	163
<i>MX 25/35 Miniature XYZ Positioners with Piezo Electric Inertial Drive</i>	164
<i>MT 25 Kinematic Mirror Tilting Stage</i>	165

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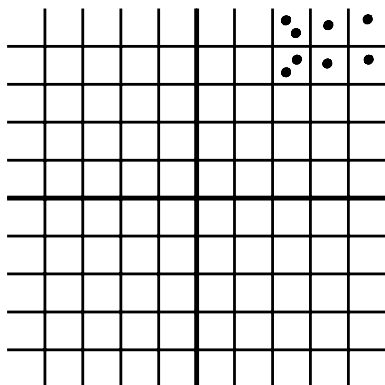
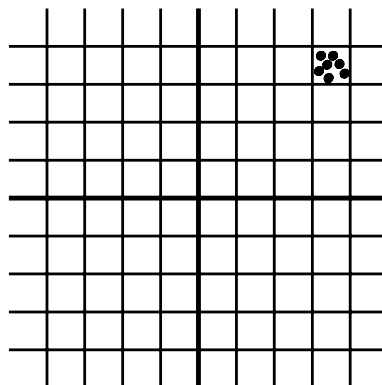
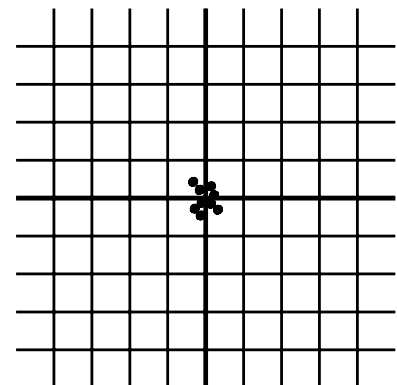
Abb. 3

Technical information

The following section explains the most important technical terms in precision positioning technology in order to make it easier for you to select the optimum positioning system for your needs.

Accuracy

The accuracy of a linear positioning system can be divided into two categories:

1. The accuracy of the path itself, referred to in technical terms as guide accuracy.
2. The linear positioning accuracy along the path.

The former is a property of the guide system (guide, ball-bearings, crossed-roller bearings, etc.) and characterizes the degree of accuracy with which motion takes place in a defined direction. The latter refers to the conversion of drive into linear motion. The limiting influences are lead-screw properties as well as encoder or other feedback mechanics.

Guide accuracy

Every object has six degrees of freedom (see Fig. 1). These are the linear motions along the three orthogonal axes X, Y and Z as well as the rotations around these axes (θ_x , θ_y and θ_z). The task of a linear guide is to restrict the motion of an object precisely to one of these axes (typically referred to as X-axis). Any deviations from rectilinear motion along the X-axis are the consequences of inaccuracy in the guide system.

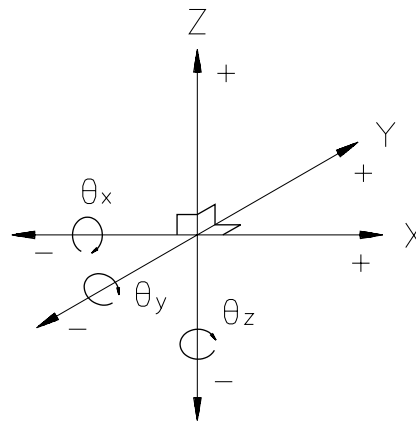
Fig. 1 Degrees of freedom

Five possible guide errors remain: motion in the X-direction, motion in the Z-direction, rotation around the X-axis (roll), rotation around the Y-axis (pitch) and rotation around the Z-axis (yaw) (see Fig. 2).

With linear positioning elements, the mainly linear motion arises from the interaction of a large number of points of support along the length of the guide. That is why any deviation of the individual guide from a straight line results both in a translational error and an angle error. As the average straightness error in the individual guide is generally extremely small, the angular errors can be neglected for most purposes. On the other hand, even small angular errors do add up over large guide lengths to considerable linearity deviations. The combination of pitch and displacement in the Z-direction yields the flatness and the combination of yaw and displacement in the Y-direction yields the straightness (see Fig. 2).

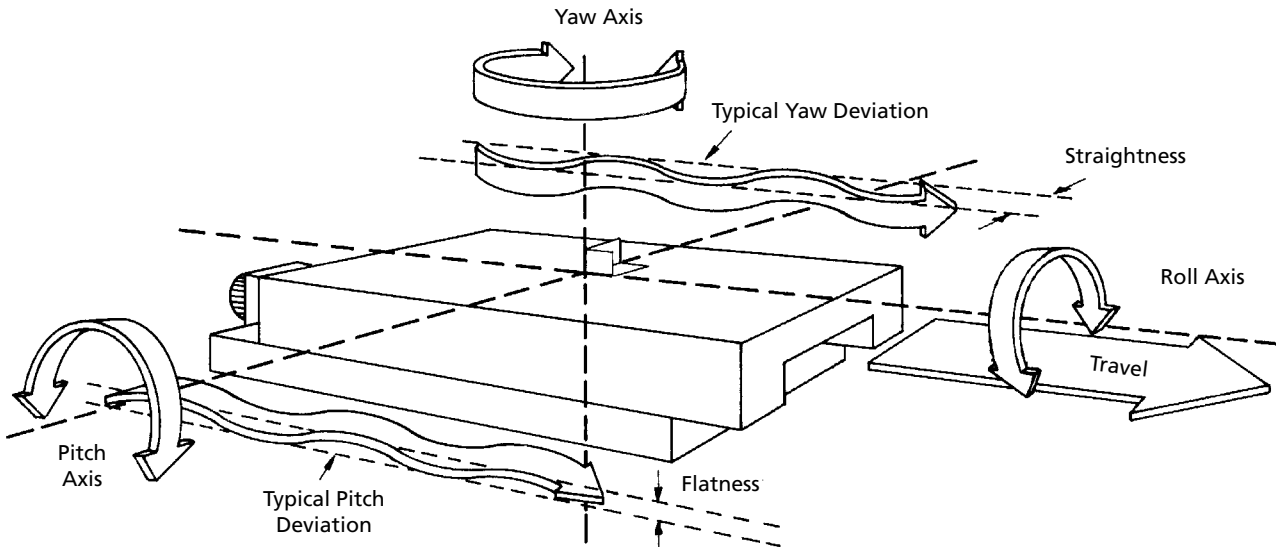
Fig. 2

Positioning Accuracy

Most linear positioning systems use leadscrews to transform the rotation of a motor into linear motion. The quality of the leadscrew determines the accuracy of the positioning system. The leadscrew nut should be free from backlash which will reveal itself upon direction reversal.

Accuracy analysis of the leadscrew error distinguishes between the cumulative error over the path and the periodic error which occurs with leadscrew rotation.

High positioning accuracy is attained using interferometric measurement, compensation data determination and online axis compensation. In practice, however, control via encoder position feedback has proved effective.

We offer two types of encoders: linear and rotary. Linear encoders eliminate a number of error sources such as leadscrew error. The accuracy of the positioning system is determined in this case by the accuracy of the encoder. Rotary encoders require considerably less space and can therefore

be more easily integrated into the positioning elements, and for that reason we use them as standard encoders.

Thermal expansion is a main source of error in high-precision positioning systems. Every degree of temperature difference to the desired temperature (usually 20 °C) results in an error of approx. $12 \cdot 10^{-5}$ due solely to thermal expansion. The thermal expansion of the positioning element itself must equally be taken into account in order to make use of its high accuracy. When using linear encoders, the thermal expansion of the encoder is the element that determines accuracy.

Reproducibility

The reproducibility of a positioning system refers to the range in which the actual position varies when a particular commanded position is to be approached. We distinguish between unidirectional reproducibility, where a position is always approached from the same direction, and bidirectional reproducibility, which refers to the approach to a position from any direction. As system-inherent inaccuracies such as nut backlash

are often not recognized with unidirectional reproducibility, bidirectional reproducibility should always be considered when there is no guarantee that the application allows a unidirectional approach to the commanded position.

A system with high reproducibility shows very little range of variation for motion to a given position, no matter which direction the approach is from.

It is important to make a clear distinction between accuracy and reproducibility. Figure 3 shows the difference between reproducibility and accuracy.

Resolution

The resolution of a positioning system is defined as the smallest position increment produced by an encoder or other feedback device. The mechanical positioning element, the drive and in some cases an electronic controller determine together the total resolution of a positioning system. In stepping motor systems, the resolution is determined for example by the leadscrew pitch.

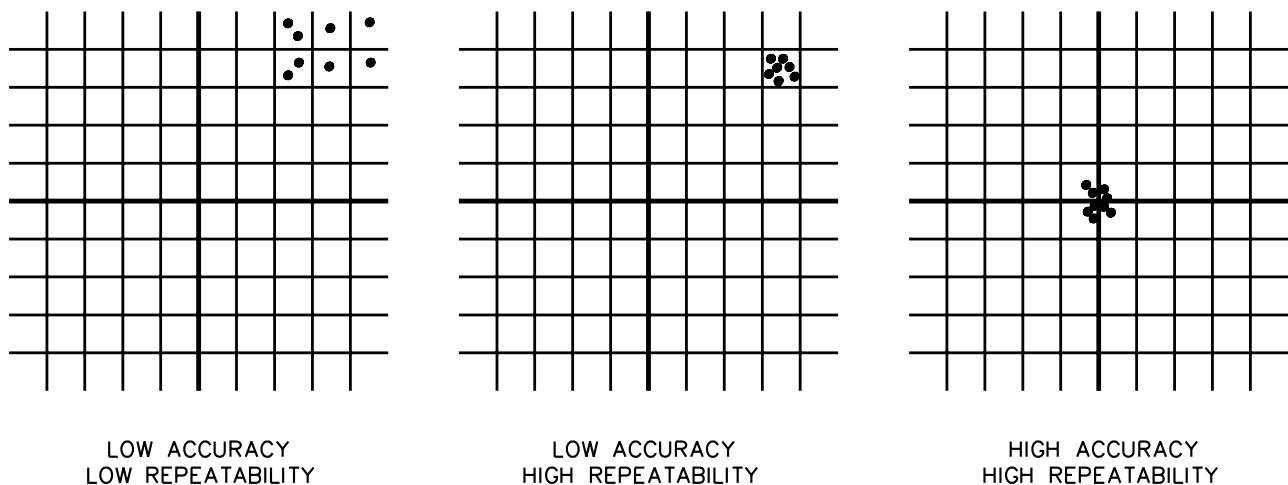
Fig. 3

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Linear Positioners TM / TMS

The linear stages TM / TMS are available in two versions, with aluminum body or stainless steel body. The sliding carriage is made of stainless steel with precision-grounded bearing surfaces for the crossed-roller bearings.


The crossed-roller bearings provide high-resolution movement and large load capacity.

These positioners can be used horizontally and vertically. They can be combined to XYZ combinations without any additional mounting brackets. Different threaded holes ensure compatibility to the LINOS construction systems.


- Compact dimensions
- High load capacity
- Large through hole
- Travel range 5 mm - 25 mm
- Crossed-roller bearings
- Micrometer with 0.25 mm pitch (TM 30-05, TM 50-16)
- Micrometer with 0.5 mm pitch (TM 90-25)
- The adapter plates allows you to combine positioners of different sizes.

Linearpositionierer TM Linearpositioner TM


Produktbezeichnung	Stellweg (mm)	Belastbarkeit im Zentrum (N)	Belastbarkeit gegen Federn (N)	Axiale Belastbarkeit gegen Mikrometerschraube (N)	Gewicht (kg)	Part No.
Item Title	Travel (mm)	Central load capacity (N)	Max. axial load against springs (N)	Max. axial load against micrometer (N)	Weight (kg)	Part No.
Linearpositionierer TM 30-05 aluminum	5	150	20	50	0.04	G403370000
Linearpositionierer TM 50-16 aluminum	16	500	20	120	0.14	G403371000
Linearpositionierer TM 90-25 aluminum	25	1000	40	200	0.48	G403372000
Linearpositionierer TMS 30-05 stainless steel	5	150	20	50	0.06	G403380000
Linearpositionierer TMS 50-16 stainless steel	16	500	20	120	0.18	G403381000
Linearpositionierer TMS 90-25 stainless steel	25	1000	40	200	0.55	G403382000
Adapter plate 90-50 aluminum						G403373000
Adapter plate 50-30 aluminum						G403374000

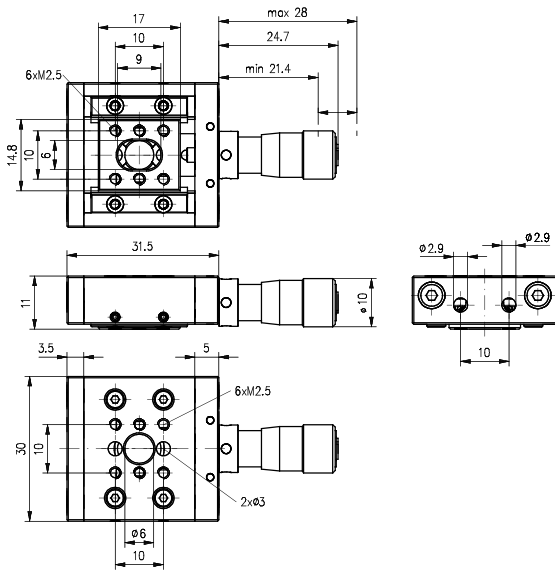Linear stage 30-05

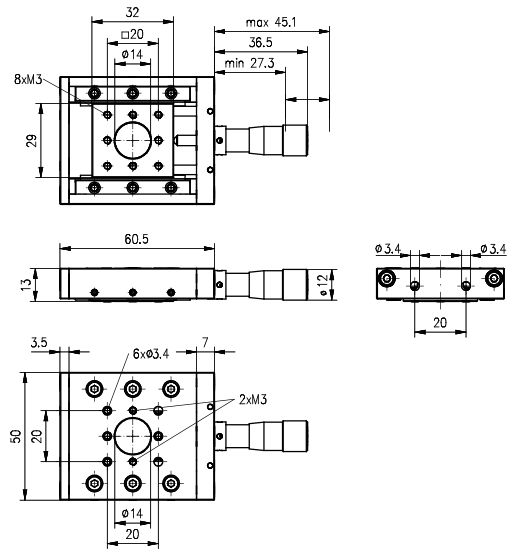
Linear stage 50-16

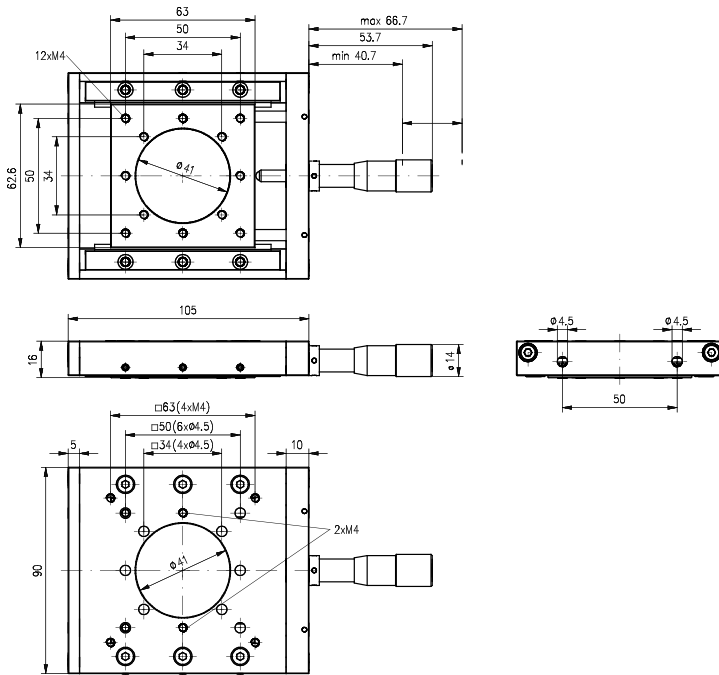
Linear stage 90-25

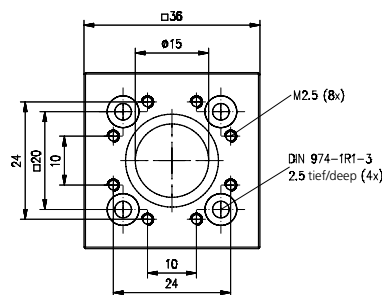
Adapter plates

Adapter plate 50-30

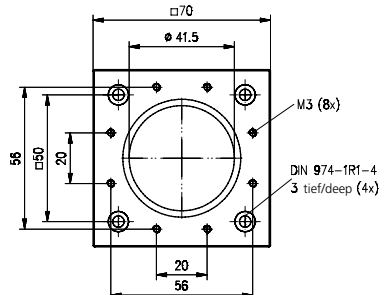Adapter plate 90-50

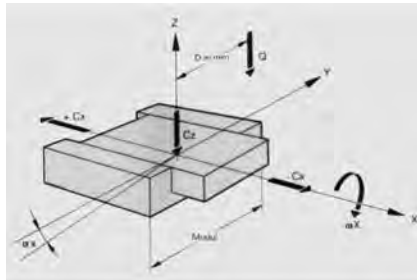Loads equivalent to the rated load Cz can be calculated using the following formula:

$C = Q (1 + D/a)$
 C = load equivalent to the central load
 Q = off-center load
 D = load distance in mm
 a = design parameter

Note that C may not assume values higher than the respective rated load Cz.

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Linear Stages TB

The TB positioner series features compact construction and high load capacity.

The ball bearing guides are bedded directly in the stainless steel base structure for exceptional guiding accuracy. Compression springs minimize backlash.

These positioners can be attached, using the threaded connectors, for use as an XY-combination. Mounting angles are available for the Z-axis.


Thanks to the large clear aperture on the positioners, they can also be used for adjustment of optical elements in the beam path.

- Travels of 5, 16, 25 and 40 mm
- Accurate movement without any play
- Ball bearing guides

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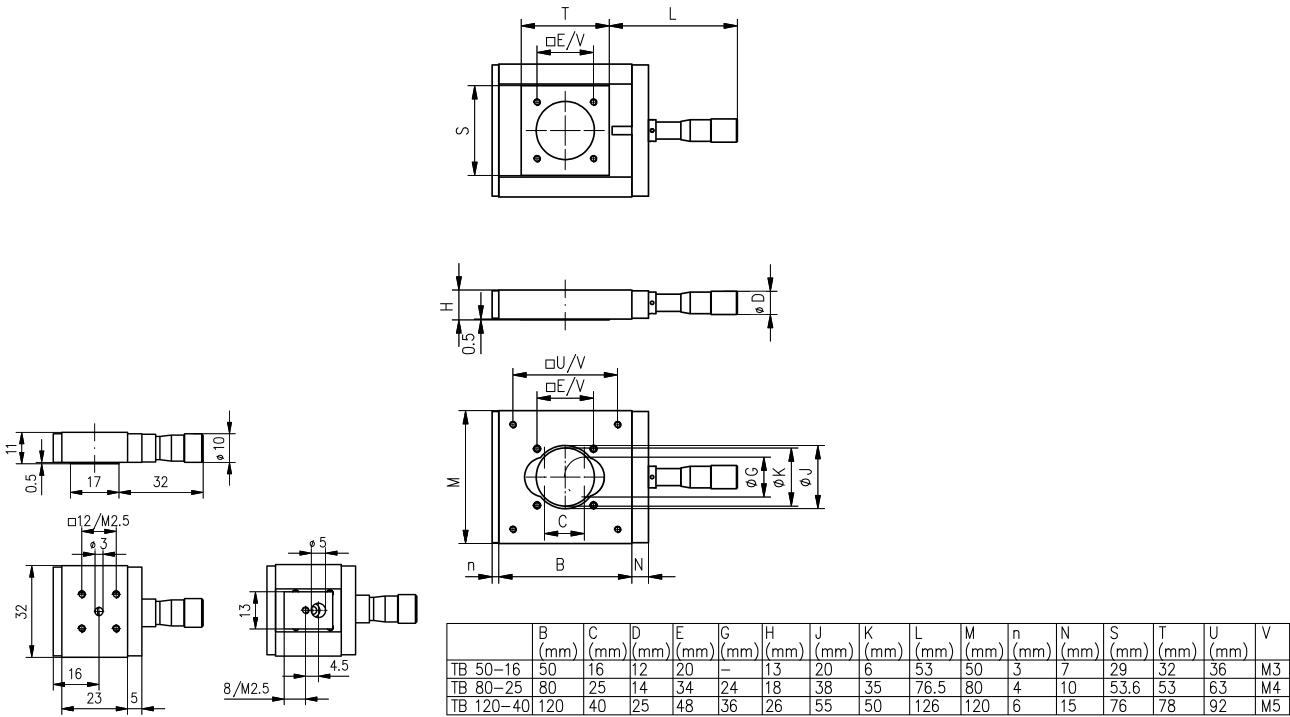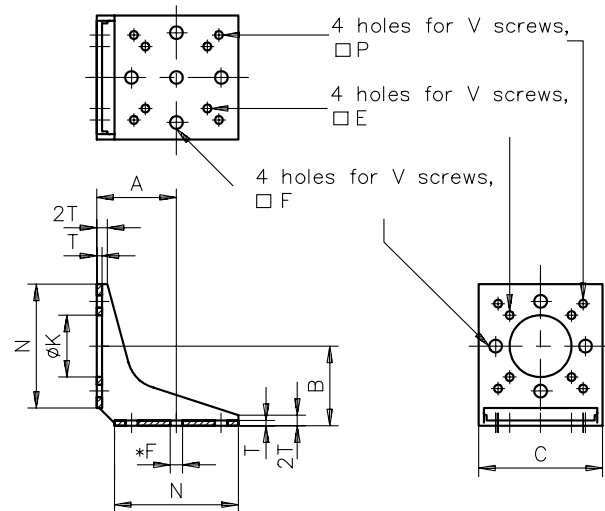
Mounting Bracket WM 50-1/80-1 for Linear Stages TB

- For external mounting of TB 50 and TB 80 linear stages

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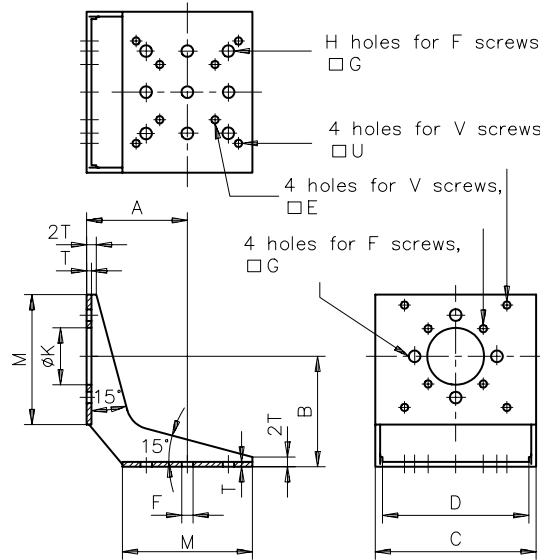
Mounting Bracket WM 50-2/80-2 for Linear Stages TB

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- For internal mounting of TB 50 and TB 80 linear stage

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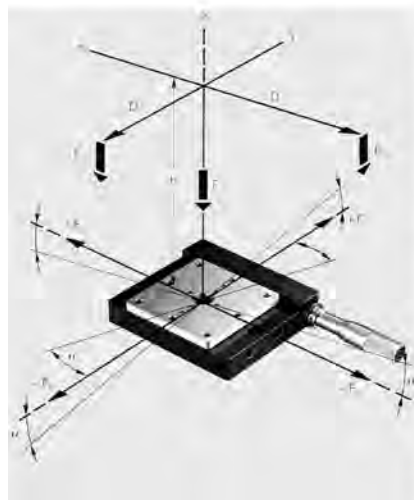
Linear Stages T 10 M/F and TS 10 M/F

The linear stages are available in four versions, central or offset drives, with micrometer screw or with fine adjustment screw.

The stages T 10 F and TS 10 F are equipped with a fine-adjustment screw with 0.25 mm pitch.

The spindle has a hexagon socket and can be driven by the provided knurled hexagon knob. If the hexagon socket is not needed, the knob can be glued in.

- High-precision ball bearings
- Accurate movement without any play
- Center (T) or lateral (TS) drive
- With micrometer (type M) or fineadjustment (type F)
- 10 mm travel
- Screw pitch 0.25 mm
- Micrometer resolution 10 µm
- Sensitivity 1 µm (type M)

Load capacity

We distinguish between two kinds of load capacity:

Rated load and maximum load.

The rated load value is an ideal weight under which all stage accuracies are maintained. The maximum load is a value which, if exceeded, could cause bearing damage.

$$F_{Dix} = \frac{F_z}{\Phi_1 + 1} \quad F_{Diy} = \frac{F_z}{\Phi_2 + 1}$$

F_z rated load, central
+ F_{sz} load on compression spring
- F_{sz} load on drive unit
± F_y lateral load

F_{Dix} off-center load, x-direction
 F_{Diy} off-center load, y-direction


α_x tilt angle, x-direction
 α_y tilt angle, y-direction

D_x load distance, x-direction
 D_y load distance, y-direction
 h height of the applied load


Φ_1 system parameter
 Φ_2 system parameter

Lineartische T 10 M/F und TS 10 M/F Linear Stages T 10 M/F and TS 10 M/F


Produktbezeichnung	Verkipfung in ax (mrad/Nm)	Verkipfung in ay (mrad/Nm)	Antrieb	Nennlast FZ (N)	Max. Belastung FZ (N)	Max. Belastung +FX (N)	Max. Belastung -FX (N)	Max. Belastung ±FY (N)
Item Title	Deflection in ax (mrad/Nm)	Deflection in ay (mrad/Nm)	Drive	Rated Load FZ (N)	Max. load Fz (N)	Max. load +FX (N)	Max. load -FX (N)	Max. load ±FY (N)
T 10 M	1.5	1.8	center micrometer	10	60	10	20	60
TS 10 M	1.5	1.8	lateral micrometer	10	60	10	20	60
TS 10 F	1.5	1.8	lateral screw	10	60	10	20	60
T 10 F	1.5	1.8	center screw	10	60	10	20	60

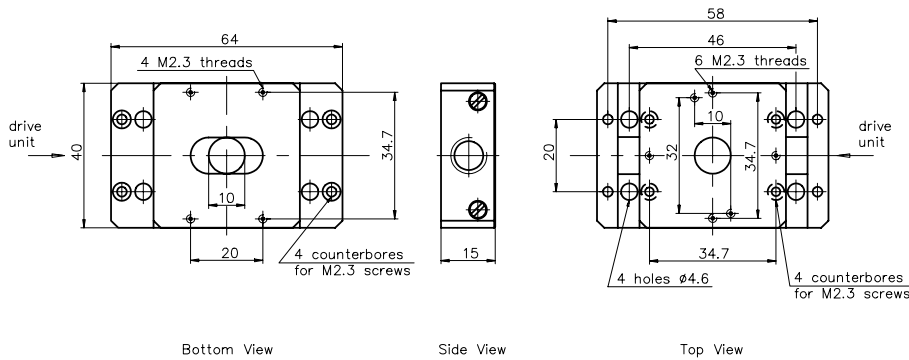T 10

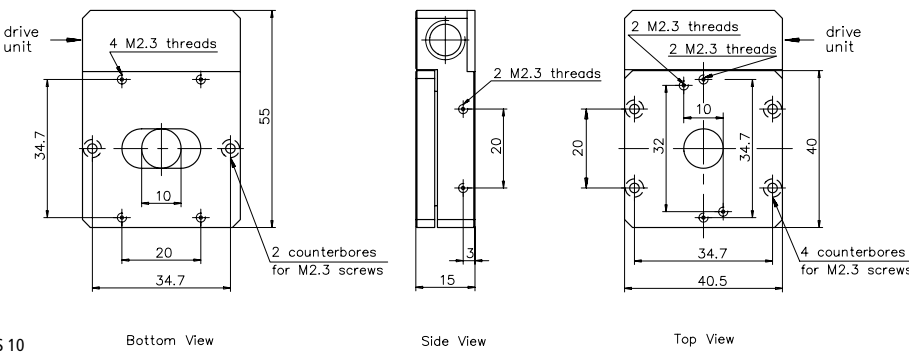TS 10

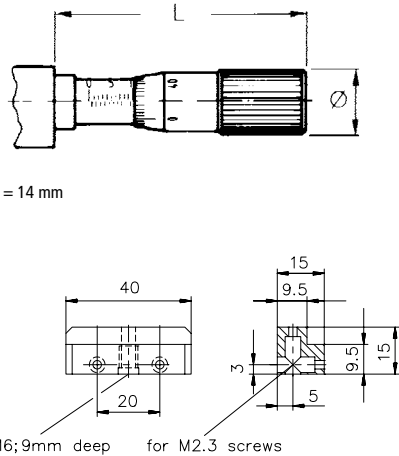
To integrate the T / TS linear stages to the Microbench.

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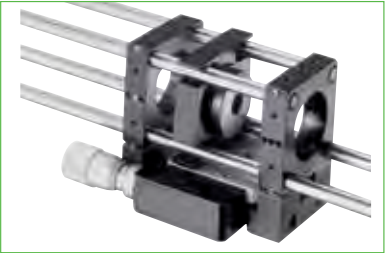The positioniers series T 10 / TS 10 are ideally suited for applications in Microbench system.

Führungs- genauigkeit unbelast. (µm)	Führungs- genauigkeit bei Nennlast F_z (µm)	Wiederholgenauigkeit unbelastet (µm)	Wiederholgenauigkeit bei Nennlast F_z (µm)	System- parameter Φ_1 (mm)	System- parameter Φ_2 (mm)	Gewicht (kg)	Part No.
Guide accuracy, no load (µm)	Guide accuracy at load F_z (µm)	Repeatability, no load (µm)	Repeatability at load F_z (µm)	System parameter Φ_1 (mm)	System parameter Φ_2 (mm)	Weight (kg)	Part No.
1	3	0.5	4	10	20	0.1	G080111000
0.5	2	0.1	1	10	20	0.1	G080211000
0.5	2			10	20	0.1	G080212000
0	3			10	20	0.1	G080113000

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

Linear Stages TL 45 / TL 65

- Preloaded ball bearing guides
- High loads
- Long life
- Micrometer resolution 10 μm
- Sensitivity 1 μm
- Mounting thread for micrometer screw M10x0.5

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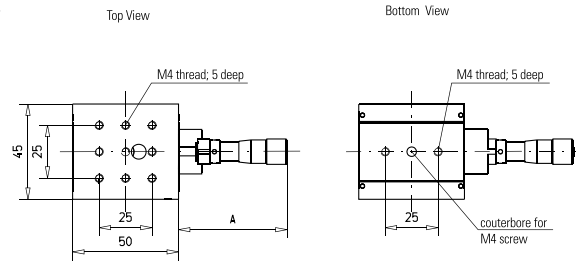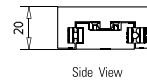
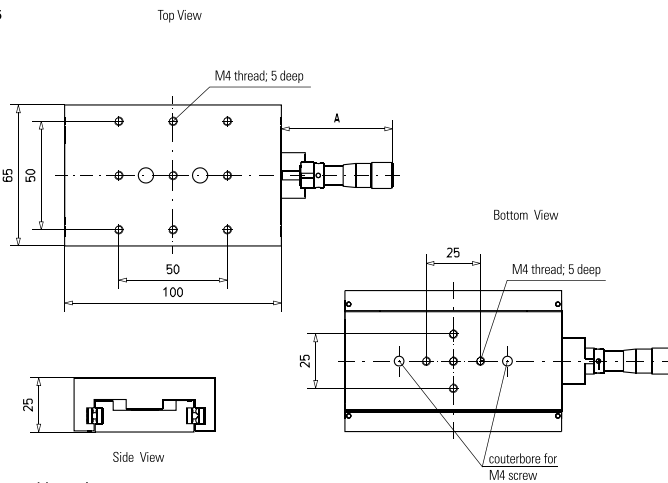
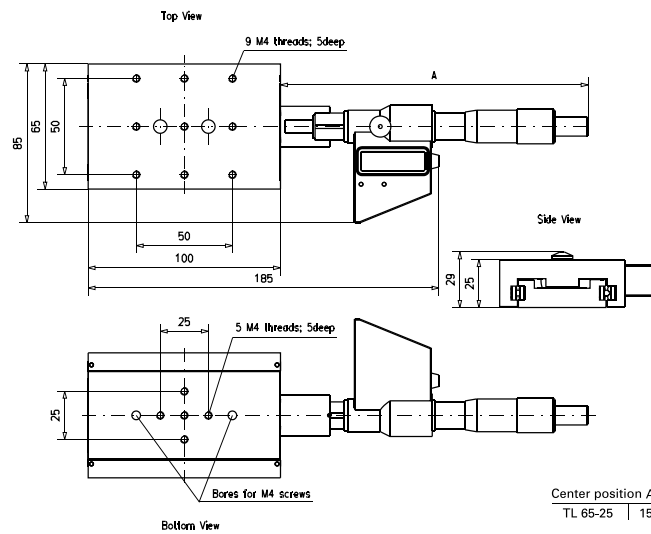
Linear Stages TL 45 / TL 65 with Digital Micrometer

- 5-digit LCD display
- Easy, correct readout
- Millimeter or inch display
- Variable zero point definition
- Absolute or incremental display
- Storage of last value
- Keeps preset values and display
- Display switched off automatically after 2 hours
- Battery operation (6 months)

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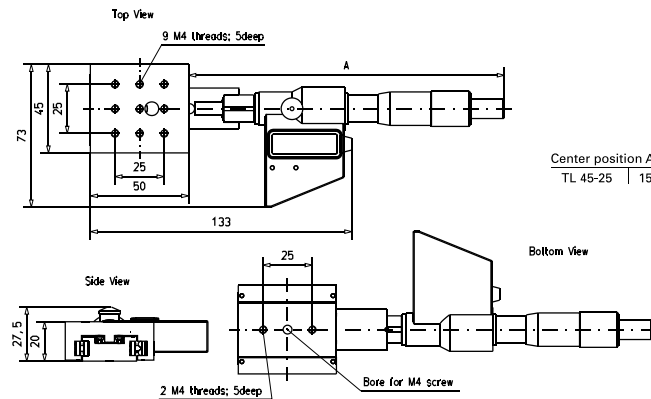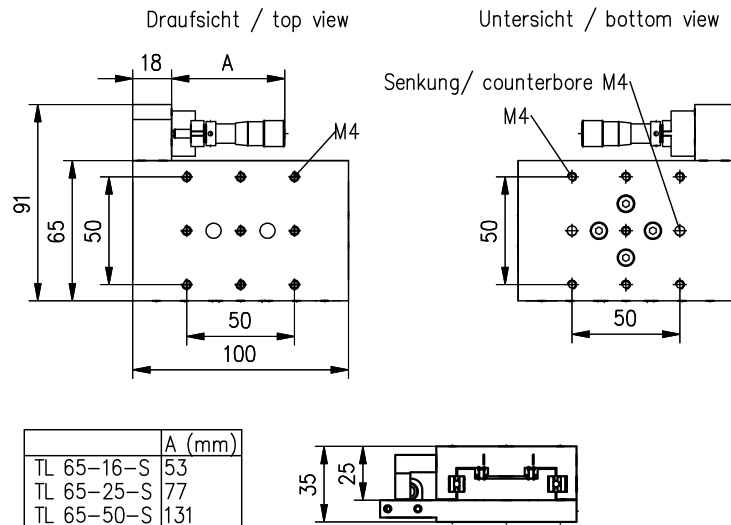
Linear Stages TL 65 with Side Drive

- Preloaded ball bearing guides
- High load capacity
- Long life
- Position is lockable
- Micrometer resolution 10 µm
- Micrometer sensitivity 1 µm

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
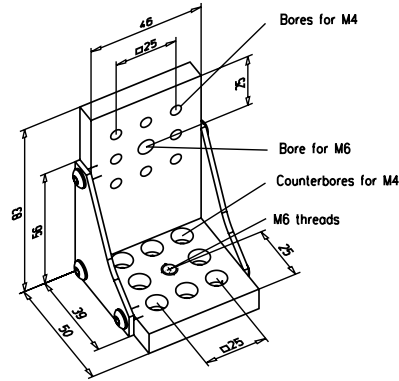
For X-Z combination of Linear stages TL-45
Compatible with Rotary stage RT 60

Mounting Components

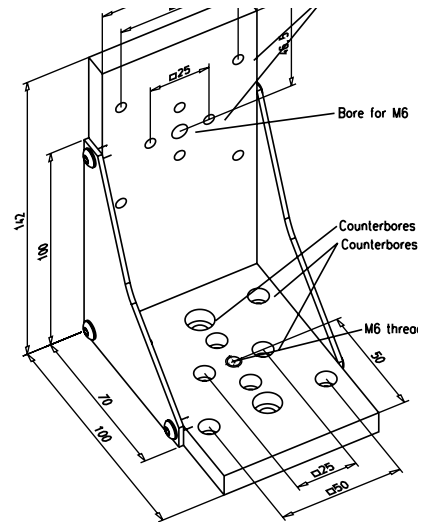
- For XZ combination of linear stages TL
- Compatible with most manual positioners

For X-Z combination of Linear stages TL-65
Compatible with stages of the XY 60/85,
RT 90/120 and HT series

Montagewinkel Z Mounting brackets

Produktbezeichnung	Part No.
Item Title	Part No.
Mounting bracket Z45	G403326000
Mounting bracket Z65	G403327000

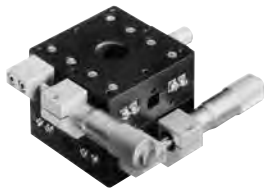XY 60-15: Extremely compact construction

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
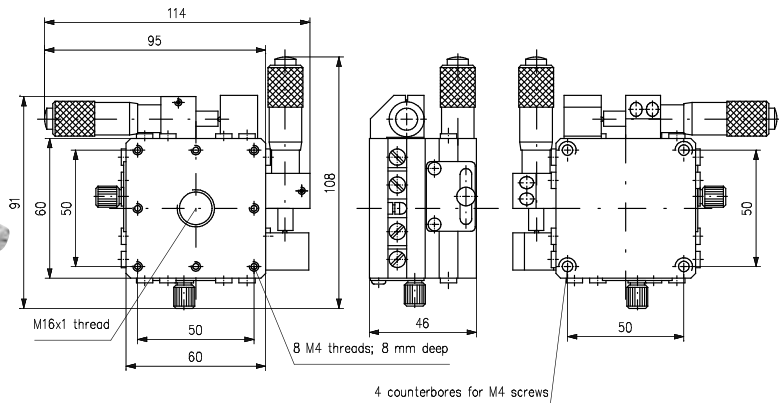
X-Y Stages XY 60 / XY 85

- Preloaded ball bearing guides
- Long life
- 10 μm scale reading
- 1 μm sensitivity

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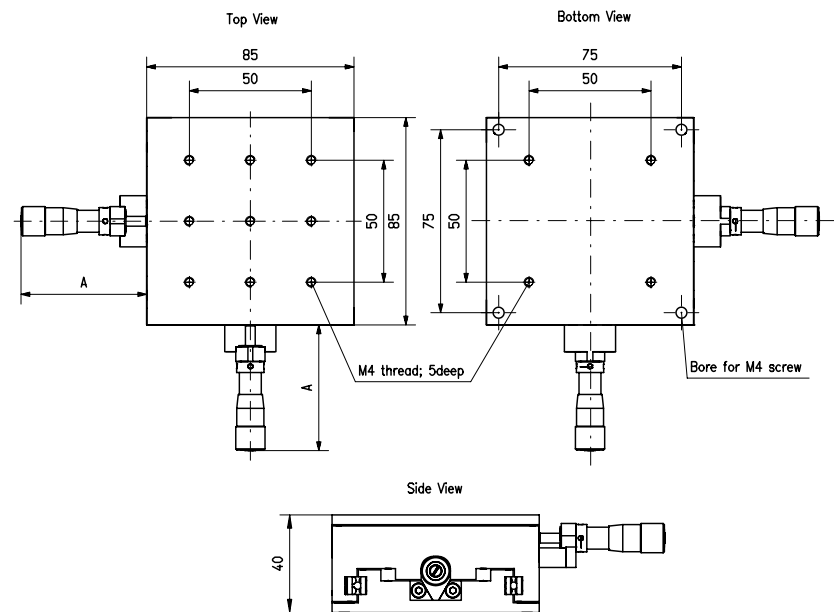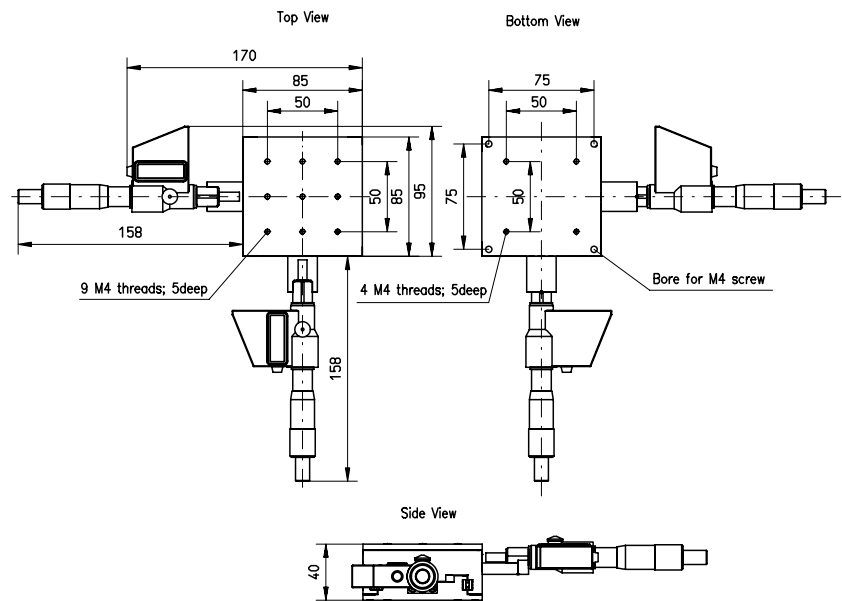
X-Y Stage XY 85-25 with Digital Micrometer

- 5-digit LCD display
- Easy, correct readout
- Millimeter or inch display
- Variable zero point definition
- Absolute or incremental display
- Storage of last value
- Keeps preset values and display
- Display switched off automatically after 2 hours
- Battery operation (6 months)

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
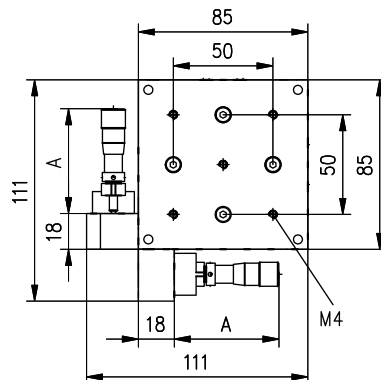
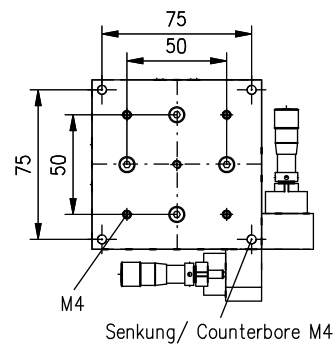
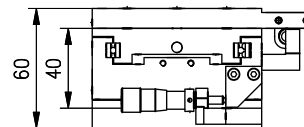
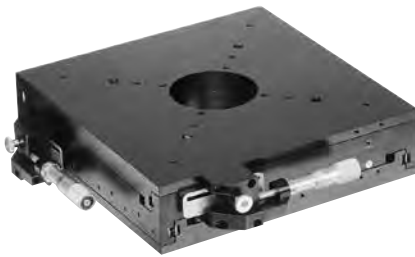
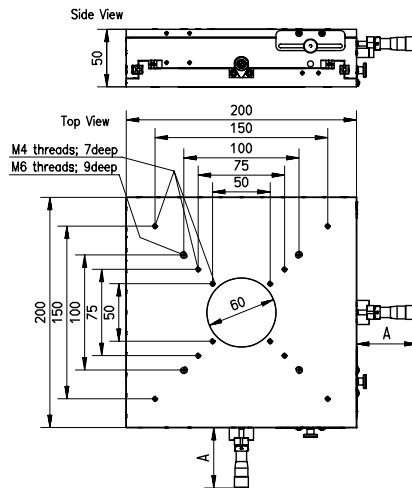
X-Y Stages XY 200


- 60 mm central hole
- Preloaded ball bearing guides
- Long life
- Position is lockable
- Micrometer resolution 10 µm
- Micrometer sensitivity 1 µm
- Side drives (-S)


Center Position A	
XY200-16	53
XY200-25	77
XY200-50	131

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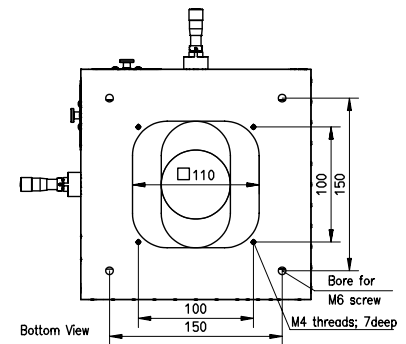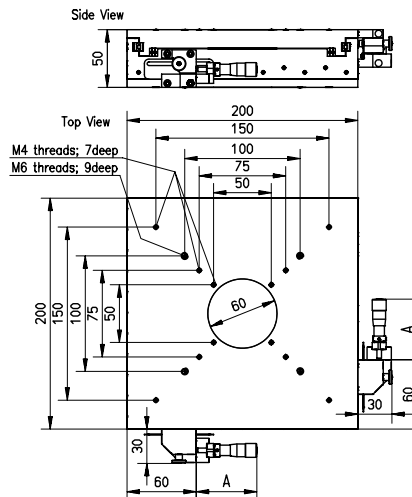
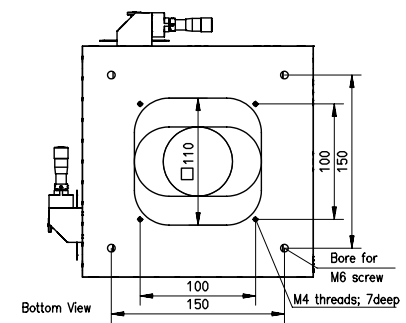
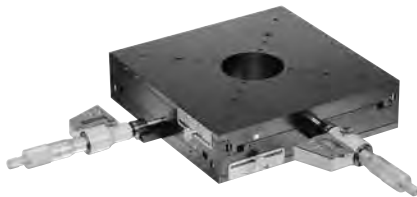
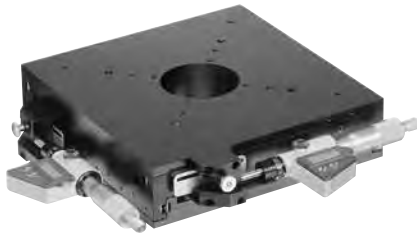
X-Y Stages XY 200 with Digital Micrometer

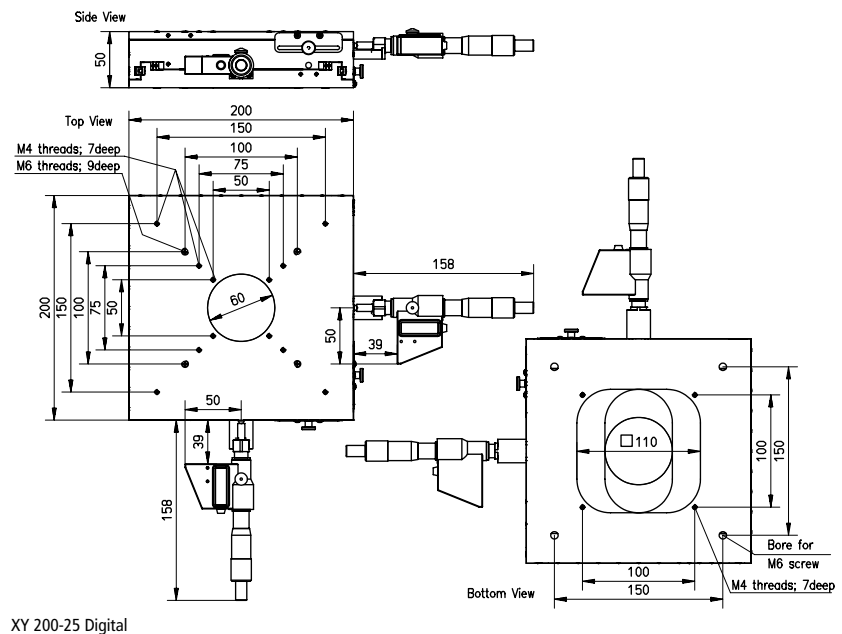- *Small overall dimensions due to sidewise mounted drives (version -S)*
- *Preloaded ball bearing guides*
- *Locking device*
- *5-digit LCD display*
- *Easy, correct readout*
- *Millimeter or inch display*
- *Variable zero point definition*
- *Absolute or incremental display*
- *Storage of last value*
- *Keeps preset values and display*
- *Display switched off automatically after 2 hours*
- *Battery operation (6 months)*

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
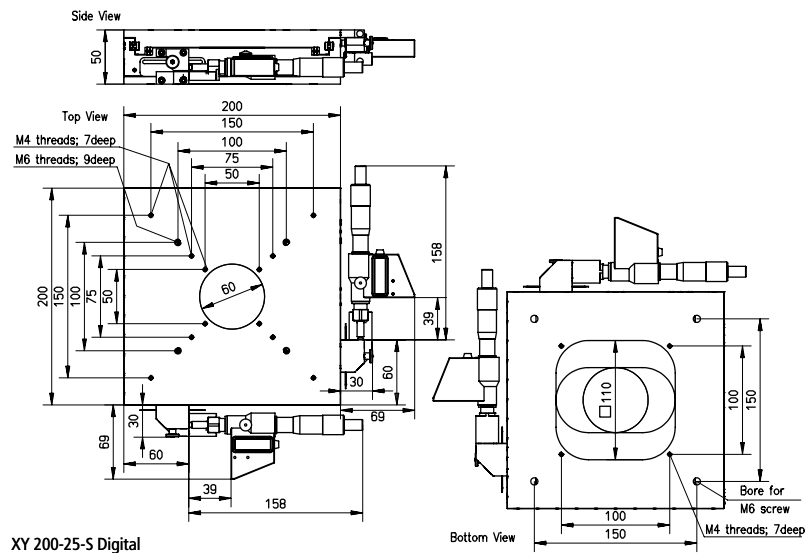
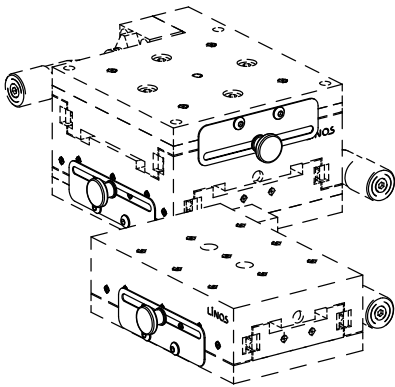
Clamping Set for Stages

- To lock position of the stages
- Suitable for linear stages TL 65 and XY Stages 85 with side drive
- Max. 50 mm travel
- 2 Sets are necessary for XY stages

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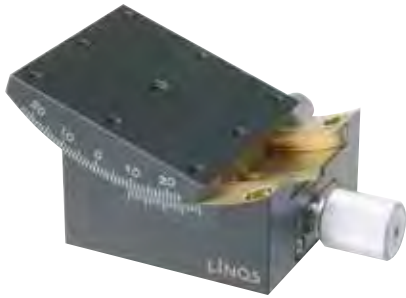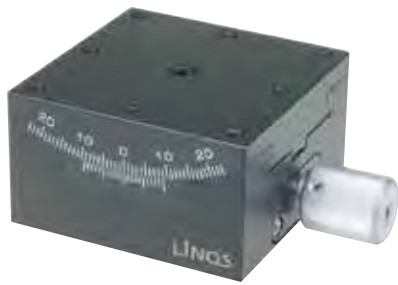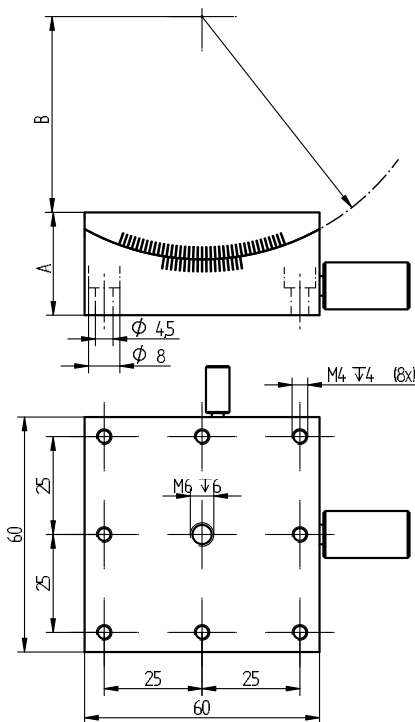
Goniometer Stages 60-60 / 60-80

Goniometer Stages enable precise angle adjustment around a fixed rotation point that is centered over the mounting surface of the positioner. You can select the radius to determine that distance to rotation point, which means LINOS Goniometer Stages can be stacked to create an adjustment range in two axes with a single rotation point.

These Goniometer Stages are made of solid brass with matte black painted surfaces and a dovetail slide for highest precision, as well as a fine adjustment screw. A vernier scale is included for reading the angle, and the angle can be fixed by tightening the adjustment screw. Compatibility with other positioners and LINOS construction systems is assured by the bore hole pattern.

- Scale with vernier
- For horizontal or vertical use
- Stage locking screw
- All adjustments are from the side
- Stages may be stacked to give true orthogonal rotation around the same external point

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
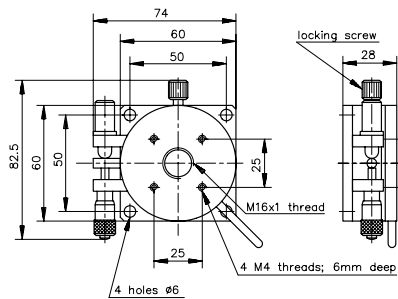
Rotary Stages RT 60 / RT 90 / RT 120

These compact Rotary Stages enable precise positioning with both coarse and fine adjustment. Coarse adjustment is performed by loosening the locking screw and then turning the rotation plate. The circumferential vernier scale lets you read off the value with a high degree of precision. Fine adjustment is performed using the laterally positioned micrometer screw. Rotary Stages can be combined with linear positioners.


- Continuous rotation
- 360°-scale reading
- Can be used horizontally or vertically
- Adjusted angle can be locked
- Laterally arranged micrometer
- Scale graduation 1°
- Parallelism 0.05 mm

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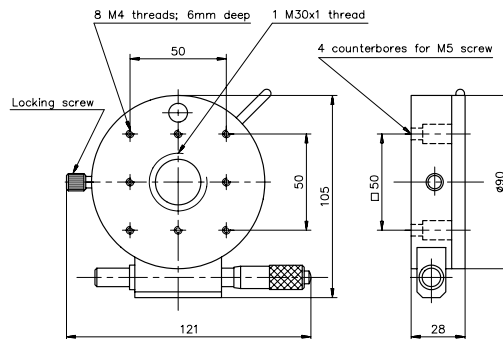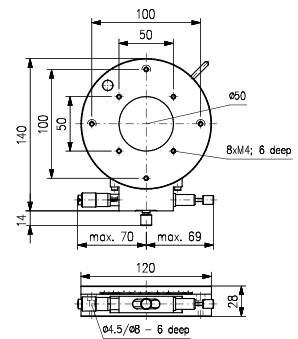
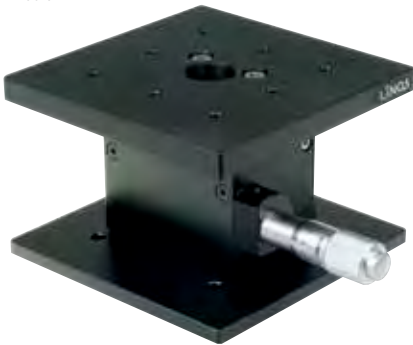
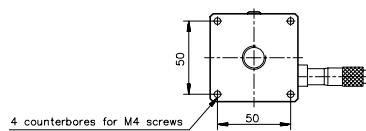
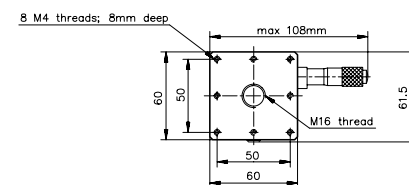
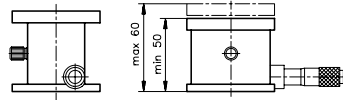
HT 60-5



Z-Axis Stages HT 60-5 / 90-5 / 120-10

These high-precision Vertical Translation Stages enable vertical positioning that is accurate to within micrometers, for optical components or similar construction elements. The micrometer screws have a readability of 10 μm and a sensitivity of 1 μm . The desired position can be fixed by tightening the clamp screw mounted on the side of the translation stage. Vertical Translation Stages are available in three sizes.

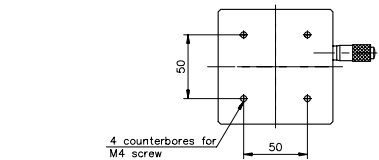
- Compact construction
- High-precision ball bearings
- Micrometer
- Locking device to secure position
- Travel ± 5 mm respectively ± 10 mm for the HT 120-10
- Sensitivity 1 μm

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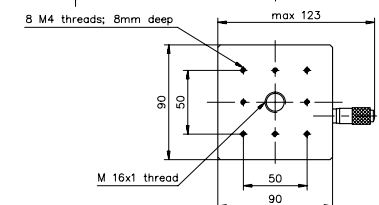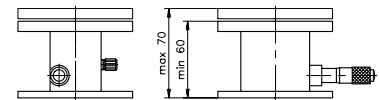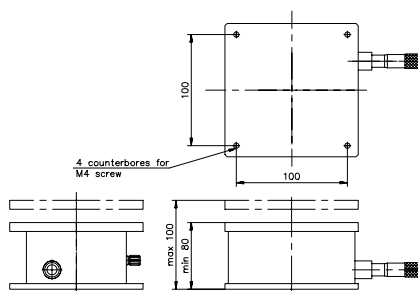
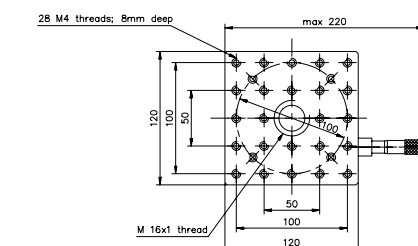
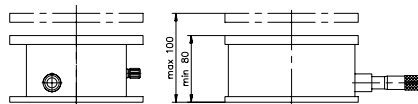
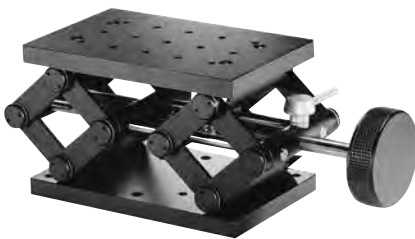
Vertical Translation Stage HT 100-40

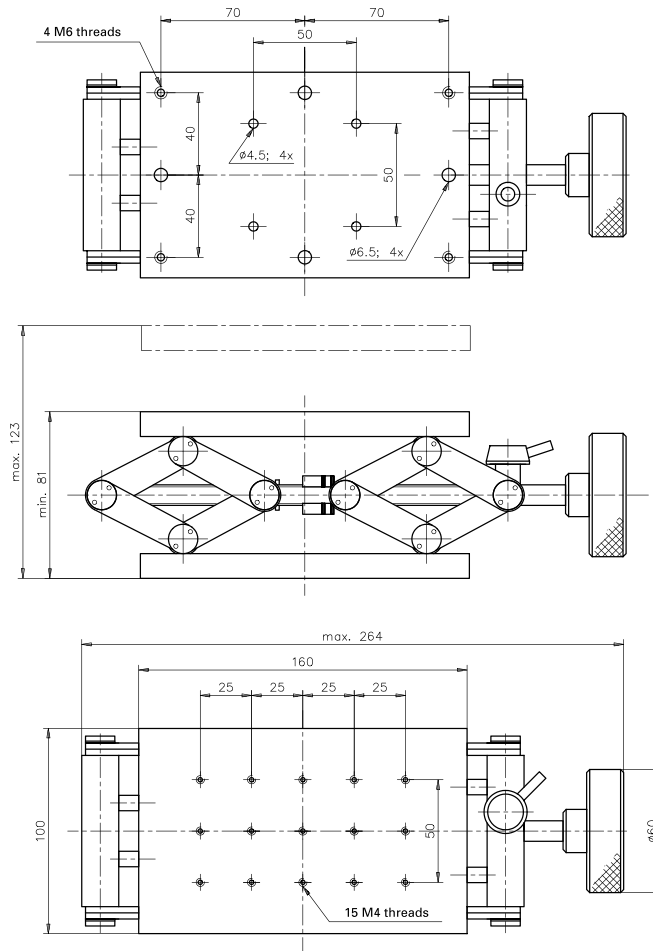- Large travel range
- Zero backlash scissors jack
- Screw rested in ball bearings
- Locking mechanism
- High load capacity

**Hebetisch
Vertical Translation Stage**

Produktbezeichnung	Stellweg (mm)	Belastbarkeit (N)	Gewicht (kg)	Part No.
<i>Item Title</i>	<i>Travel (mm)</i>	<i>Load Capacity (N)</i>	<i>Weight (kg)</i>	<i>Part No.</i>
Vertical Translation Stage HT 100-40	40	100	1.6	G402184000

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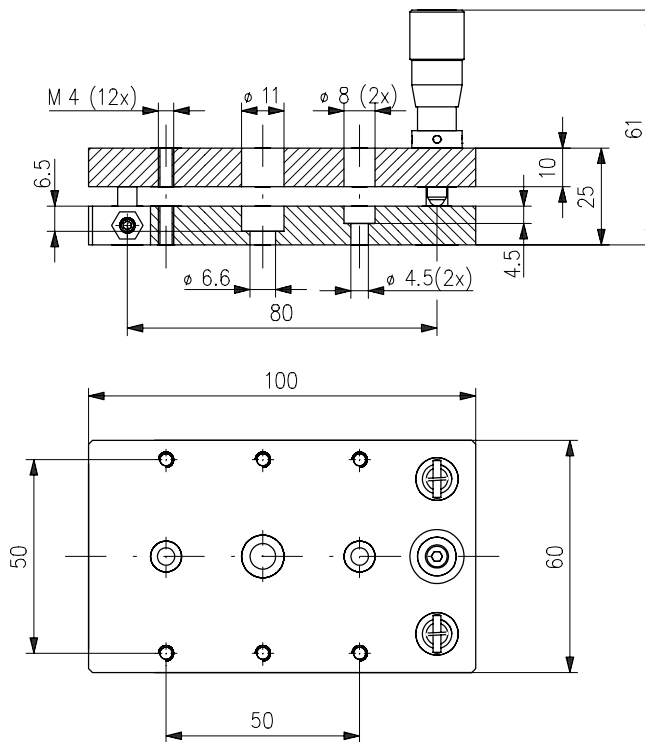
Tilting Stage NT 60-3

The Tilting Stage enables precise tilting, even for heavy components. It gives you an adjustment sensitivity of 3" with a range of $\pm 3^\circ$.

This positioner is made of high-strength aluminum with a matte black anodized surface for resiliency and scratch resistance.

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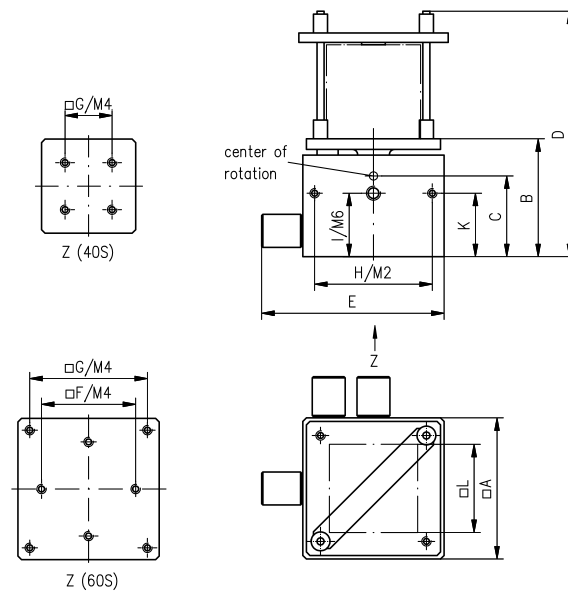
Prism Stages 40 S / 60 S

These Prism Stages let you perform adjustment along three axes, the ideal adjustment technique for optical parts such as prisms or beam-splitter cubes. The optics are gently fastened to the mounting platform with a bracket; no adhesive is necessary. The optics can be adjusted within in three of degrees freedom around a common center of rotation. The Prism Stages come with fine adjustment screws as standard equipment. These screws can be exchanged by micrometer screws for model 60 S, upon request. This lets you read the position you set.

The height of the optical axis of the prism stage 40 S is 40 mm. This ensures the compatibility with Microbench.

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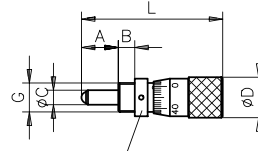
Locking collar
(Tool: G06 1098 000 Hook wrench)

Micrometers G

- Fits to Holder for micrometer or Rod holder G06 1229 000 (both chapter Microbench)
- Locking collar to adjust zero position
- 10 µm resolution

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Precision Adjustment Screws G, bushing with connection thread

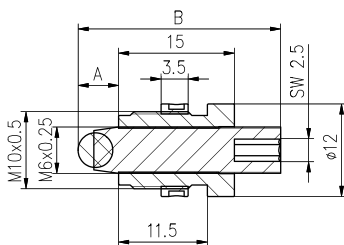
The fineadjustment screw G consists of a stainless steel spindle in a brass bushing. That ensures a smooth run. The pitch of the thread is 0.25 mm. The screws are available with flat tip or ball tip.

The spindle has a hexagon socket and can be driven by the provided knurled hexagon knob. If the hexagon socket is not needed, the knob can be glued in.

The fineadjustment screws can be mounted to many LINOS elements by using the M10x0.5 outer thread with locking collar.

Please order the hook wrench separately.

- 0.25 mm pitch leadscrew
- hexagon socket
- knurled knob included

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
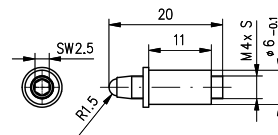
Fine-adjustment screws M4

Fine-Adjustment Screws, Flat Bushing

The fine-adjustment screws consists of a stainless steel spindle in a brass bushing. That ensures a smooth run. The pitch of the thread is 0.15 or 0.25 mm. The screws are equipped with ball tip. The spindle has a hexagon socket.

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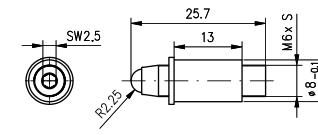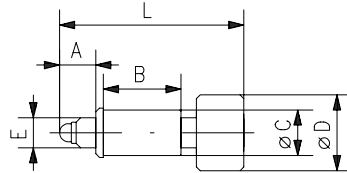
Precision Lead Screws Lees

- Lees™ technology
- Smooth adjustment
- Very high resolution

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Datasheets for all Piezo Positioners can be found on the respective product pages at www.qioptiq-shop.com below the "Docs + Drawings" button.

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CN 30 Controller for Piezo Electric Inertial Motors

- *Special design for piezo driven step motor*
- *Works with batteries or external power supply*
- *Operates up to 3 axes*
- *Applications without position sensor or closed loop*

Controller for Piezo Electric Inertial Motors

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Piezo-Controller CN 30	G403513000

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CF 30 Controller for Piezo Electric Inertial Motors

- *Special design for piezo driven step motor*
- *Works with USB-voltage or external power supply*
- *Operates up to 3 axes simultaneous*
- *Applications without position sensor or closed loop*

Controller for Piezo Electric Inertial Motors

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Piezo-Controller CF 30	G403518000

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Example: MS 30 and MX 25 mounted onto flat rail FLS 40

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
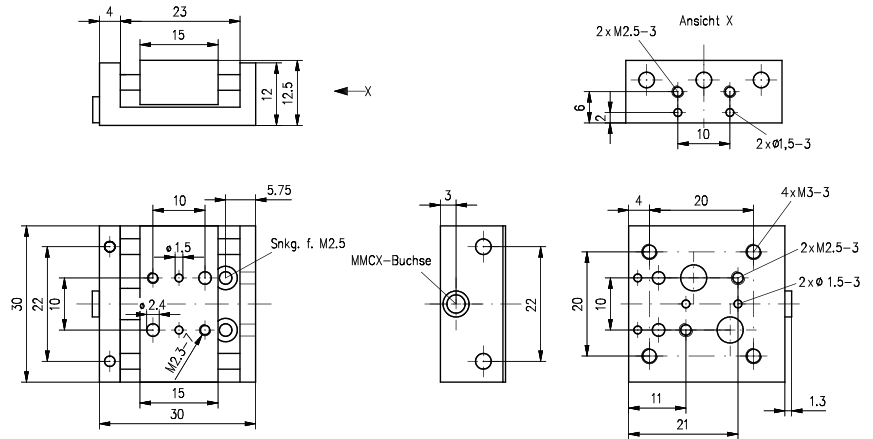
MS 30 Miniature Translation Stage

- XY or XYZ combinations
- CNC-machined aluminium body
- Precision ball bearing guides
- No limit switches necessary
- Customized designs possible
- Driven by hand-held (CN 30) or USB controller (CF 30)

- Piezo driven step motor with low hysteresis
- Holds reached position without current
- Step width about 15 nm (depends on controller)
- Velocity up to 1.5 mm / s (depends on controller)

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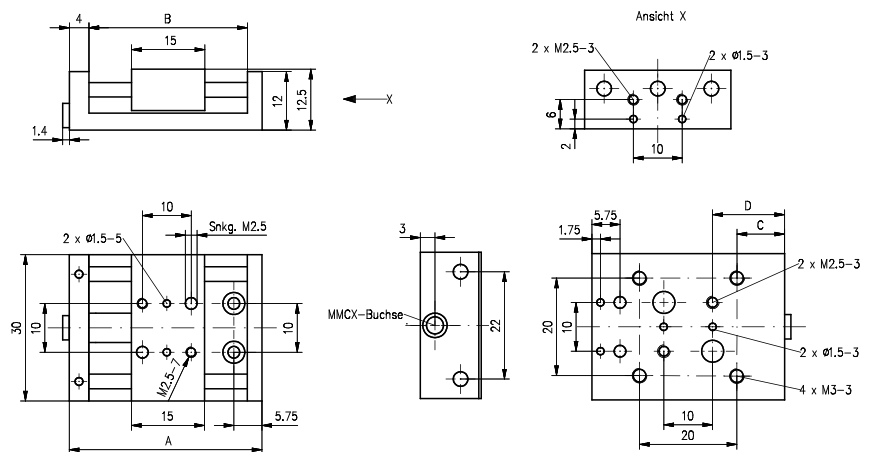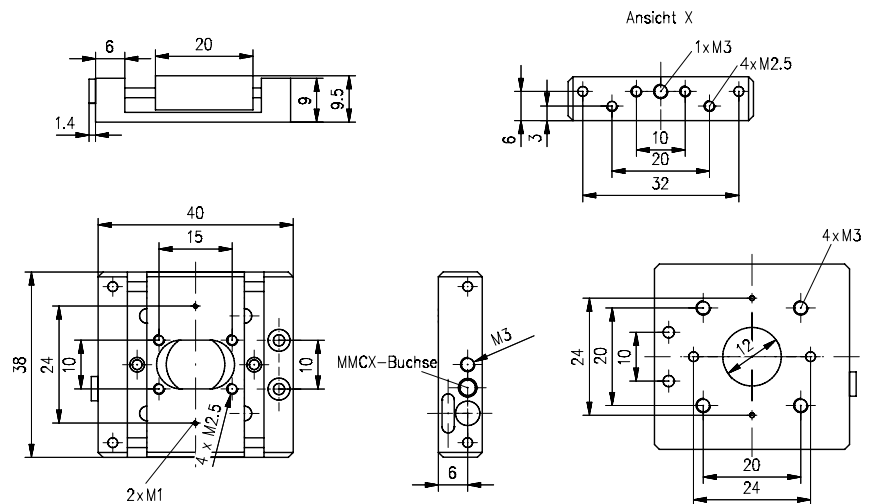
MS 38 Miniature Translation Stages with Free Opening

- Fits to optical bench systems (FLS 40)
- XY or XYZ combinations
- CNC-machined aluminium body
- Precision ball bearing guides
- No limit switches necessary
- Customized designs possible
- Driven by hand-held (CN 30) or USB controller (CF 30)
- Piezo driven step motor with low hysteresis
- Holds reached position without current
- Step width about 15 nm (depends on controller)
- Velocity up to 1.5 mm / s (depends on controller)
- Travel 8 mm
- Free opening of 10 mm (on whole travel)

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

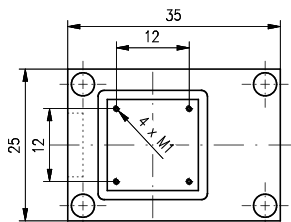
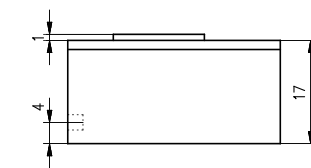
MX 25/35 Miniature XYZ Positioners with Piezo Electric Inertial Drive





- CNC-machined aluminium body
- No limit switches necessary
- Customized designs possible
- Driven by hand-held (CN 30) or USB controller (CF 30)
- Piezo driven step motor with low hysteresis
- Holds position without power
- Up to 10 mm travel in xyz
- Step width about 15 nm (depends on controller)
- Velocity up to 1.5 mm/s (depends on controller)

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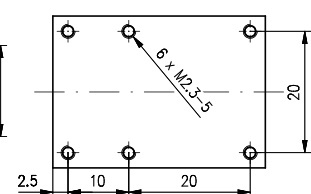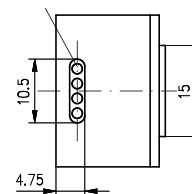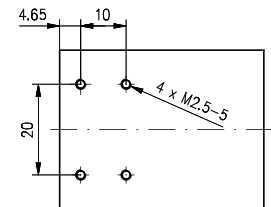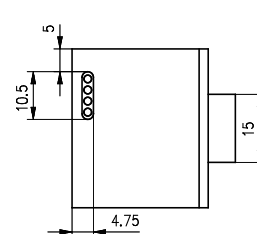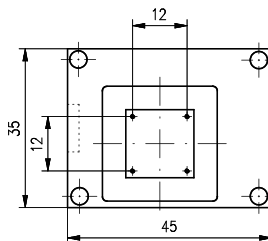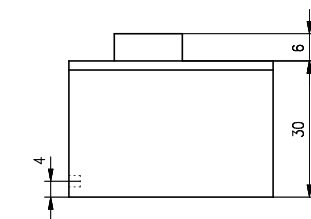The outer diameter of 25 mm enables use of the MT 25 as a motorized prism table and mirror mount in the Microbench.

MT 25 as mirror positioner in microbench setup

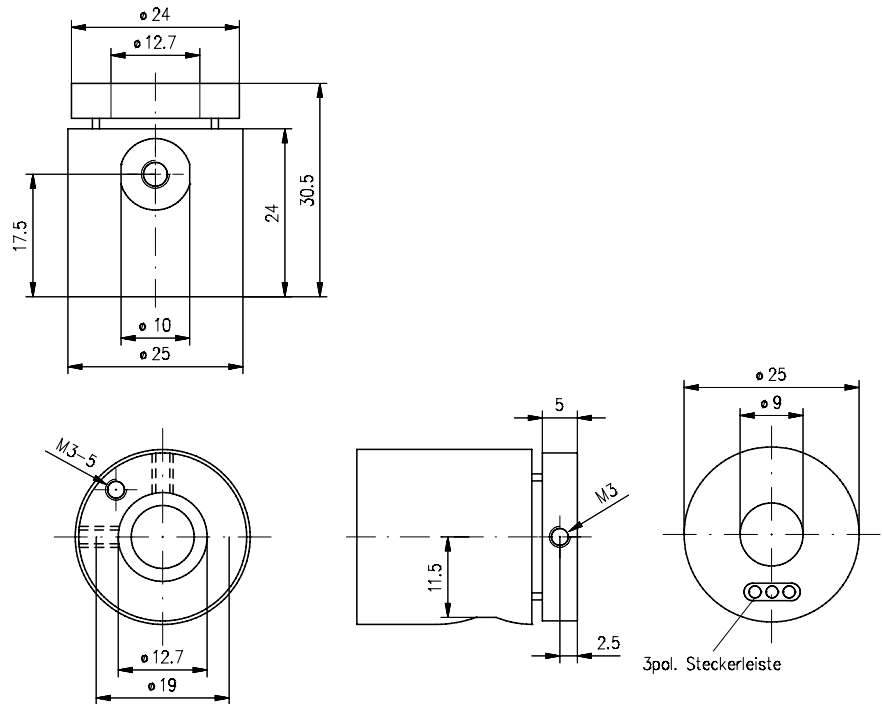
MT 25 as prism stage in microbench cube 25

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MT 25 Kinematic Mirror Tilting Stage



- Tilt in two axis
- Also usable as prism stage
- No limit switches necessary
- Customized designs possible
- Driven by hand-held (CN 30) or USB controller (CF 30)
- Piezo driven step motor with low hysteresis
- Holds reached position without current
- Optical height 11.5 mm (on rod)
- Angular adjustment 6 deg (±3 deg)
- For mirror with dia. 12.7 mm
- Free opening of 9 mm
- Solid state hinges
- Step width about 14 µrad
- Velocity up to 25 mrad/s (> 1 deg/s) (depends on controller)

MT 25 Kinematische Spiegeljustierung Kinematical Mirror Tilting Stage with piezo electric inertial drive

Produktbezeichnung	Part No.
Item Title	Part No.
Mirror Tilt MT 25	G403508000

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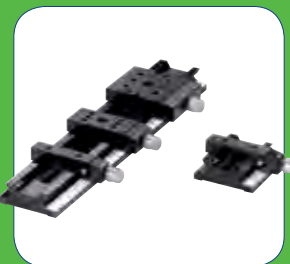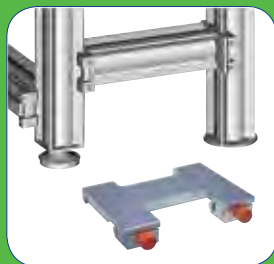Our quality criteria:

- *Precise bearing surfaces*
- *Excellent straightness; low torsion*
- *Stable design*
- *Mounting independent of existing hole patterns*

Rail Systems

LINOS flat rail systems can be used as a low-profile micro-optical bench and beam steering system. The bearing surfaces on the low-warpage aluminum alloy are precision-milled and treated with abrasion-resistant black anodizing, making these rail systems an ideal basis for linear and areal structures.

We offer three sizes, all compatible with the LINOS construction systems, to cover every application: The Flat Rail System FLS 40 is compatible with the LINOS Microbench and the FLS 95 with the LINOS mounts, posts and holders.

In addition to these flat rails, LINOS manufactures the X95 profile system especially for three-dimensional structures. All LINOS rails have slotted holes for fastening them to optical tables with M6 or 1/4" bore hole patterns. The sophisticated design, consisting of a precise rail and compatible carriers with spring-loaded clamping, ensures an extremely precise positioning of components in the structure.

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Special Features:

Carriers with springs enable high-precision repositioning of components on the rail.

Ideal Areas:



Flat optical benches in research and education; beam-steering systems; flexible assembly of components within instruments.

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
Content

Rail System 40	
Rail FLS 40	170
Carrier FLR 40	172
Rail System 65	
Flat Rails 65	174
Carrier 65	175
Profile- and Rail System X95	
X 95 Profile System	176
Carriers FLR 95	178
FLS 95 Rail System	179
Carriers X 95	180
Carrier Positioner X 95 - 30	181
X 95 Clamping Carrier	181
Cross Carrier X 95	182
Carrier X 95 - MB	182
X 95 Base Plate	183
Clamps SP 4 and SP 4-10	183
X 95 Locking Caster	184
X 95 Levelling Foot	184
X 95 End Plate	185
X 95 Corner Connector	185
X 95 Cover Plate	186
X 95 Single Leg	186
X 95 Double Leg	187
X 95 Hinged Connector	187

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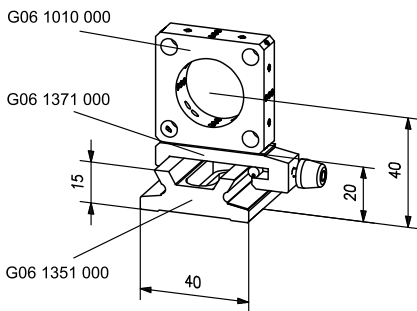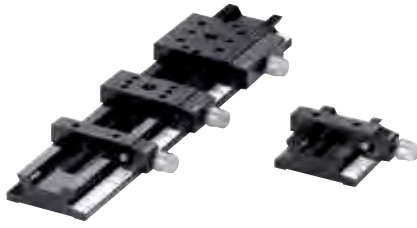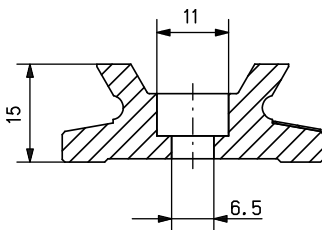Rail FLS 40

The black anodized profiled rail FLS 40 features high linearity and low torsion. Thanks to additional, precise machining, all rail and support surfaces of the extruded profiled rails are exceptionally even. Slots in the profiled rails enable components to be assembled in a choice of metric or inch standard tapped hole patterns, and the rail positions to be easily shifted, independently of the tapped pattern. A ruler makes it easy to locate the components to ensure the reproducibility of their positioning.

The FLS 40 Rail system has a high degree of levelness due to the precise processing of all guide and bearing surfaces. The rail profile has elongated holes so that it can be attached to work surfaces with metric or inch standard threads as well as independent of the pattern.

All FLS 40 rails have slotted holes for fastening them to optical tables with M6 or 1/4" bore hole patterns. The sophisticated design, consisting of a precise rail and compatible carriers with spring-loaded clamping, ensures an extremely precise positioning of components in the structure. The FLS 40 rail system is available in various lengths from 40 mm to 1,000 mm

- Flat, micro-optical bench
- High stability
- Precise guide rail surfaces
- Including ruler
- Slots allow rails to be shifted to any position independently of the tapped hole pattern
- Abrasion-resistant black anodized surface
- Material: aluminum alloy, black anodized
- Mass: 1.0 kg/m
- Straightness: 0.3 mm/m
- Torsion: 0.2°/m

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
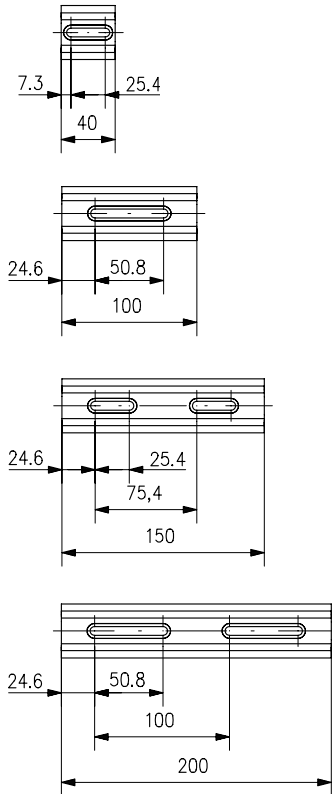
For higher flexibility you can directly combine Rail system FLS 40 and numerous Systems and components like Microbench, Laser, Positioners. Please find further information about application examples, guidelines, frequently asked questions and more in our Brochure "Microbench, Bench and Rail systems". You will find the brochure for download at each Microbench product under "Docs + Drawings".

Positions of slotted holes of rail lengths 40 up to 200 mm


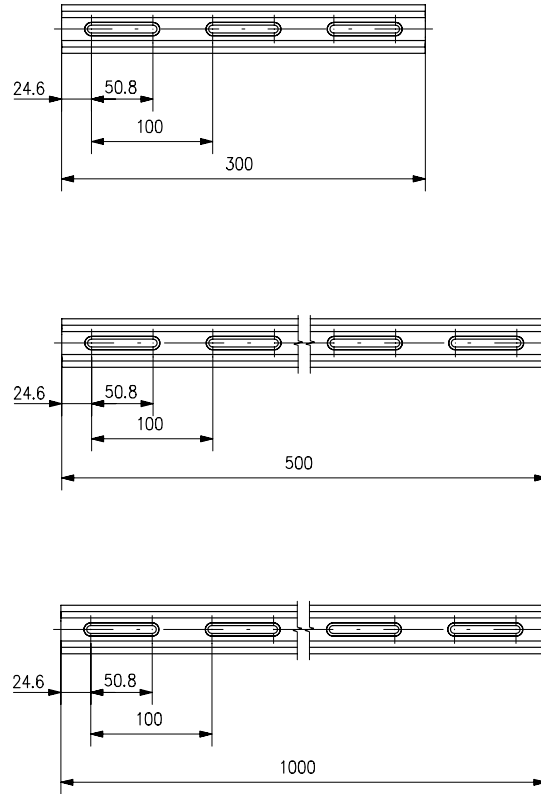
Positions of slotted holes of rail lengths 300 up to 1000 mm

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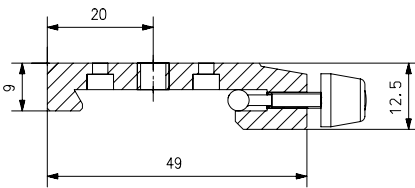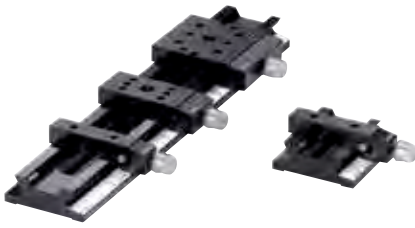Carrier FLR 40

The carriers FLR 40 are delivered in different sizes designed to connect the FLS 40 rail system to the most diverse of our system components. The carriers can be secured in place using a spring-loaded thumbscrew. When the thumbscrew is loosened, the carriers maintain their position because of the spring tension. However, they feature smooth gliding action for easy positioning. As a result, they prevent jolts when their position is shifted.

A 20 mm tapped hole pattern enables a number of components of the LINOS Microbench system to be mounted, such as mounting plates and adjustment elements. Moreover, cubes 25 and 30 of our Microbench can be mounted on the special pattern available on the FLR40-40 carrier, making it easy to assemble 3-D set-ups. A central M6 thread allows rods, posts and columns to be mounted.

- Spring-loaded clamps for smooth gliding
- Thumbscrew for securing in place
- 20 mm slot pattern for mounting Microbench components
- Optical axis height of 40 mm for use of Microbench components
- Central mounting of posts, columns and holders using M6 mounting thread
- Black anodized


statement

"We highly appreciated the increased flexibility due to various threads. Furthermore, the knurled-head screws are secured against losing."


A customer

Reiter FLR 40 Carrier FLR 40


Produktbezeichnung	Breite (mm)	Part No.
Item Title	Width (mm)	Part No.
FLR 40-10	10	G061371000
FLR 40-20	20	G061372000
FLR 40-34	34	G061373000
FLR 40-40	40	G061374000
FLR 40-80	80	G061378000

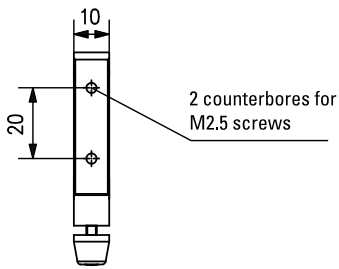
FLR 40-10 for assembling a Microbench mounting plate or Microbench positioning systems

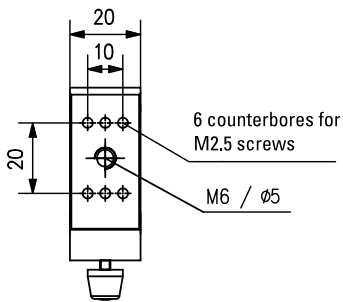
FLR 40-20 for assembling Microbench positioning systems and mounting plates as well as rods (diameter 14 mm) Nanobench base plate N, and Nanobench holder N

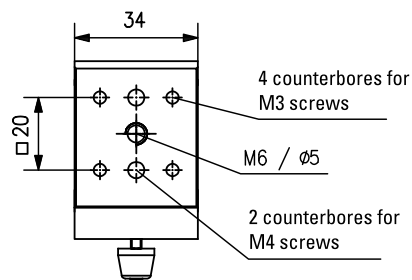FLR 40-34 for assembling Microbench positioning systems and mounting plates as well as rods (diameter 14 mm), posts 14, Nanobench holder N and mirror holder for diameter 50 and 63 mm

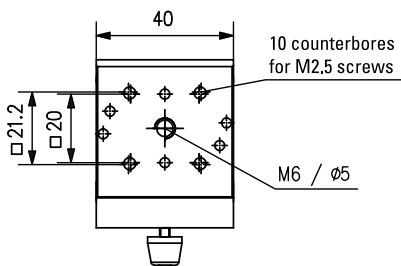FLR 40-40 for assembling Microbench cubes 25 and 30, Microbench positioning systems and mounting plates as well as rods (diameter 14 mm), Nanobench base plate N, and Nanobench holder N 20

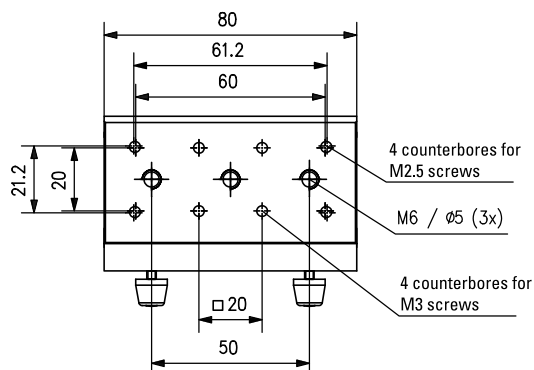
FLR 40-80 for assembling linear stages TL 45 and TL 65, Microbench cubes 25 and 30, Microbench positioning systems and mounting plates as well as rods (14 mm diameter), post 14, Nanobench base plate N, and Nanobench holder N 20

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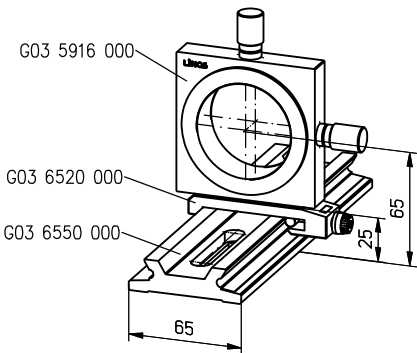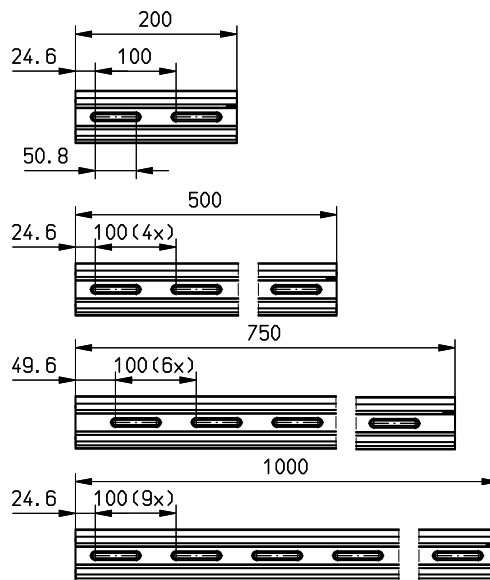
Flat Rails 65

The flat, extruded rail features excellent straightness properties and low torsion. Bearing surfaces are precision processed for smooth movement of the carrier. The rail features a matte black anodized, abrasion-resistant surface.

Slotted holes make it easy to fasten the rail to optical tables independent of existing hole patterns, using either M6 or 1/4" screws.

The FLS 65 rail is available in lengths from 200 mm to 1000 mm.

- Material: aluminum alloy, black anodized
- Mass: 1,8 kg/m
- Straightness: 0,6 mm/m
- Torsion: < 0,2°/m

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


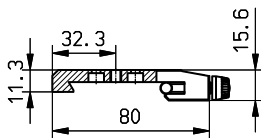
Carrier 65

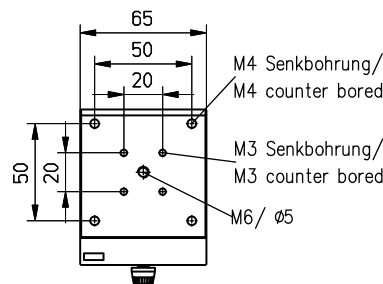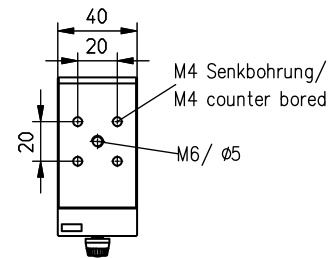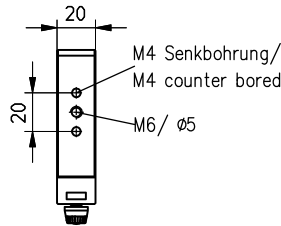The carriers are available in three widths and feature a universal bore hole pattern for attaching optical mounts. The spring-loaded clamp ensures smooth movement of the carrier along the rail, and a knurled screw fixes the carrier in position. Like the rails, the carriers also feature matte black, abrasion-resistant anodizing. Together with the rail, the surface of the carrier is 25 mm above the base.

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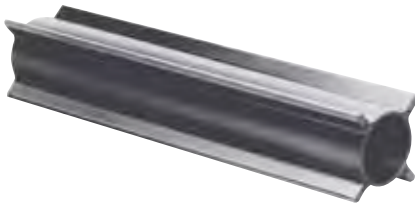
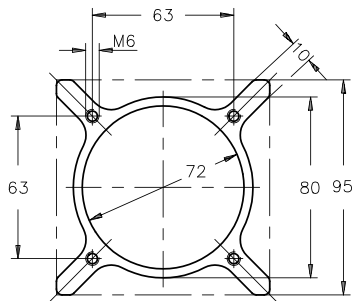
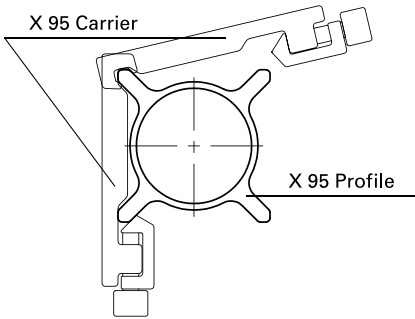
X 95 Profile System

The X95 profile system includes both profiles and carriers, and is a versatile construction system for optics and precision mechanics. It is particularly well suited as an optical bench, or for table and frame structures.

An X95 Profile is a hollow, cylindrical aluminum profile with four lateral ridges which reinforce the system and also hold the carriers. These ridges increase the torque of inertia and torsional strain considerably. Four threaded bore holes on the profile end planes provide fastening options for connectors or caps. The X95 profiles are available in many standard lengths. Custom lengths of up to 3 m can be supplied on request.

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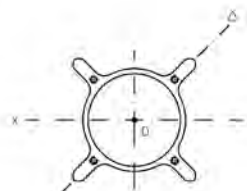
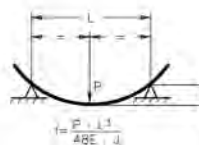
- For assembling tables or framework structures
- Hollow extrusions with cruciform arrangement of longitudinal ribs
- Allows assembling of three-dimensional structures
- Smooth-gliding carriers equipped with spring-loaded clamps
- Easy assembling
- Mount on vibration-isolation platforms/columns
- Anodized Aluminum
- Moments of inertia $J_x / J_{\Delta} \sim 220 \times 10^4 \text{ mm}^4$
- Torsional stress $J_0 \sim 1.2 \text{ kg cm}^2$ per cm length
- Young's modulus $E \sim 70 \text{ kN/mm}^2$
- Tolerance Warpage 0.5 mm/m
- Tolerance Twist 0.6 mm/m

X 95 Profile strength/dimensional-tolerance data

Moments of inertia $J_x / J_{\Delta} \sim 220 \times 10^4 \text{ mm}^4$
 Torsional stress $J_0 \sim 1.2 \text{ kg cm}^2$ per cm length
 Young's modulus $E \sim 70 \text{ kN/mm}^2$

Tolerances
 Warpage: 0.5 mm/m
 Twist: 0.6 mm/m²

f Deflection
 P Transverse load
 L Unsupported length
 J Moment of inertia
 E Young's modulus

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Please find further information and application examples in our Brochure "LINOS X 95 System". You will find the brochure on www.qioptiq-shop.com for download at all X 95 products under "Docs + Drawings".

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The stable base for vertically mounting X 95 profiles is our Mounting base, G024343000. The base is equipped with levelling feet, additionally the feet allow a height adjustment up to 20 mm. Please find the Mounting base described in chapter "Mounts and Posts", section "Posts, Columns and Rods".

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
Carriers FLR 95

- Simple carriers for flat rail system FLS 95 and profile system X 95.
 - Black anodized aluminium.
 - Fastening of the carriers on the profile by a clamping screw with plastic cone point.
 - Without spring loaded clamps, no moving without backlash possible.
- Carriers width 28 mm or 80 mm
 - Carriers with spring loaded clamps see carriers X 95.

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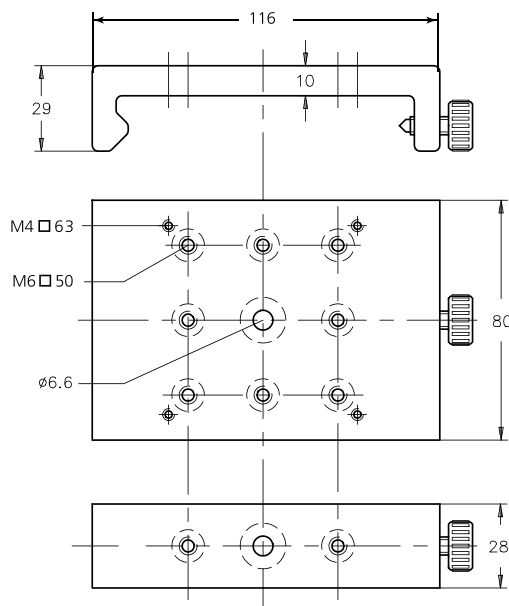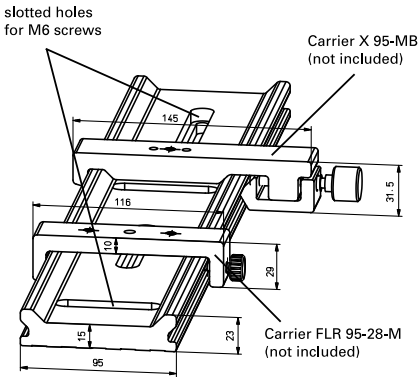
FLS 95 Rail System

The Rail System FLS 95 is ideal for use as a low-profile optical bench, particularly for setups on optical tables. Its slotted holes make it independent of existing hole patterns, and fit both M6 and 1/4" screws. The profile corresponds to the X95 system and can use the same carriers. The rail surface is treated with abrasion-resistant black anodizing and the bearing surfaces are milled to ensure the highest guiding accuracy.

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
- Low-profile optical bench
- Ideal for mounting components or instruments
- Slots in rails allow for readily shifting rail positions
- Extremely rigid
- Rails accept X 95-system components

- Aluminiumprofil, schwarz eloxiert

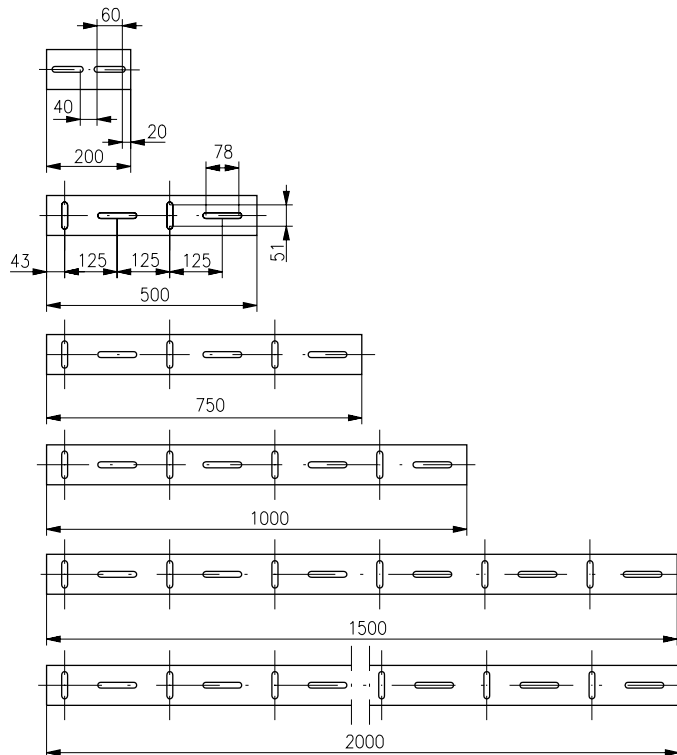
- Black anodized profiled aluminum rails

Locations and positions of slots for M6 mounting screws

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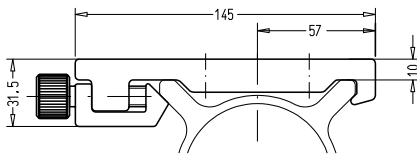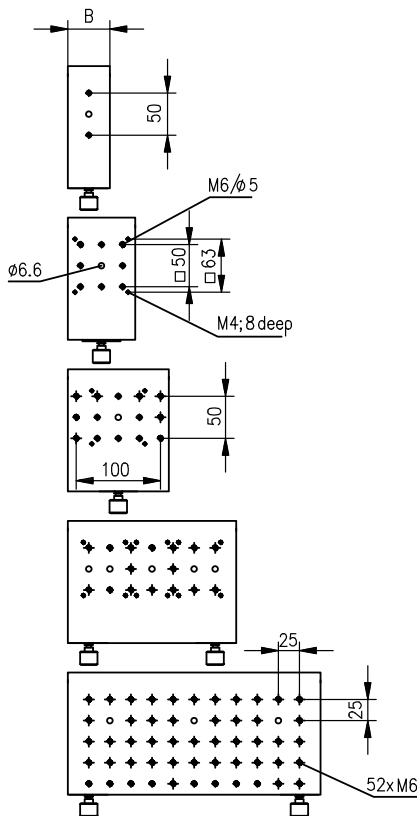
Carriers X 95

- Smooth gliding due to spring-loaded clamps
- Teflon-coated surfaces
- Anvil/clamping screw locking mechanisms
- Easy handling
- With or without hole pattern

- \varnothing 6.6 mm center bore for posts or similar devices
- M6 threaded hole for top mounting, 50 mm hole pattern
- Bottom mounting with counterbored holes for M4 screws, 50 mm hole pattern
- M4 mounting threads, 63 mm hole pattern, for mounting plate MP 80 and mounting bases

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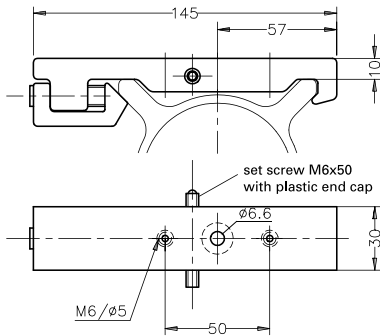
Please view chapter "Mounts, Posts and Holders" for a complete range of compatible components, such as:

- Optic and Object holder
- Plate holder and Mounting plates
- Posts and Columns

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
Carrier Positioner X 95 - 30

- For positioning X 95 Carriers (travel: up to 20 mm)
- For mounting X 95 Double leg, G02 6303 000 to X 95 Profile
- \varnothing 6.6 mm center bore for posts or similar devices
- M6 threaded hole for top mounting, 50 mm hole pattern
- Bottom mounting with counterbored holes for M4 screws, 50 mm hole pattern

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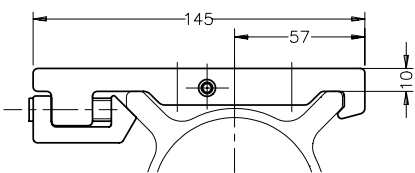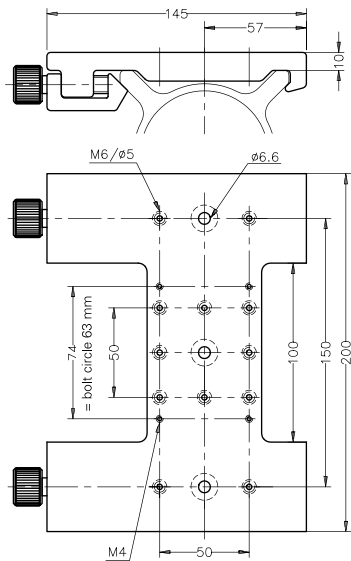
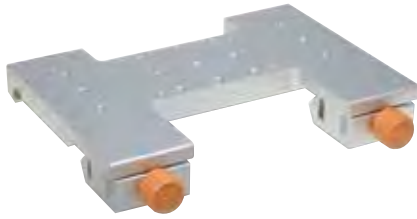
CAD files of all mechanical components can be found on the respective product sheets at www.qioptiq-shop.com below the button "Docs and Drawings" .

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
Cross Carrier X 95

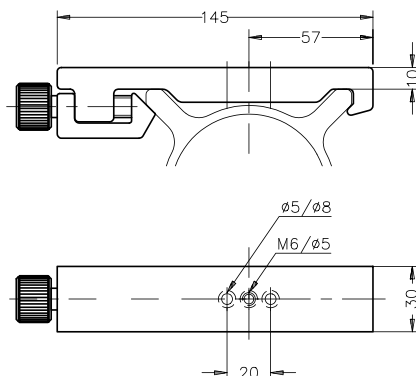- For mounting components at the intersection of two optical axes such as beam steering mirrors or beamsplitters
- Teflon-coated guide surfaces
- Anvil / clamping screw locking mechanisms
- Two anvil clamps for rigid mounting
- Equipped with spring-loaded clamps for smooth gliding actions
- $\varnothing 6.6$ mm center bore for posts or similar devices
- M6 threaded hole for top mounting, 50 mm hole pattern
- Bottom mounting with counterbored holes for M4 screws, 50 mm hole pattern
- M4 mounting threads, 63 mm hole pattern, for mounting plate MP 80

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Carrier X 95 - MB

- For bottom mounting of Microbench mounts
- Teflon-coated guide surfaces
- Anvil/clamping screw locking mechanisms
- Equipped with spring-loaded clamps for smooth gliding action

- For bottom mounting of Microbench mounts
- Mounting of posts, etc. via central M6 thread
- With counterbored holes for M4 screws

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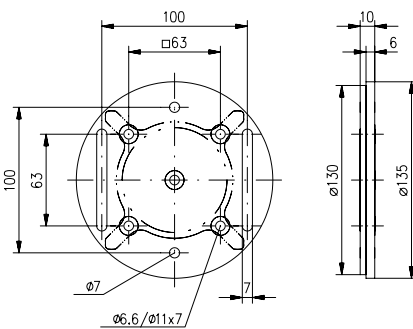
X 95 Base Plate

For perpendicular mounting of profiles on optical tables.

- Easy positioning with slotted holes
- Supplied complete with mounting screws
- Easy clamping with Clamps SP 4
- Colorless anodized aluminum

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Clamps SP 4 and SP 4-10


For clamping items on FLR or X 95 carriers. For clamping base plates or angle brackets having suitable tapped-hole patterns M4 square 63 mm.

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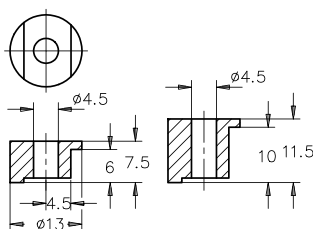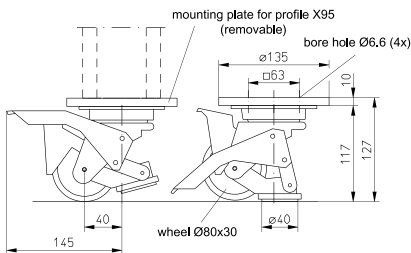
X 95 Locking Caster

For use on mobile X 95 profile structures

- Caster with integrated, stable foot
- Easy to use locking lever
- Load capacity of up to 2 kN
- Mounts on X 95 corner connectors or ends of X 95 profiles
- Supplied complete with mounting screws

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
X 95 Levelling Foot

- Removable dowel allows mounting of breadboards
- Laminated vibration isolator plate diameter 100 mm, 15 mm thick
- Maximum load capacity of 3.5 kN

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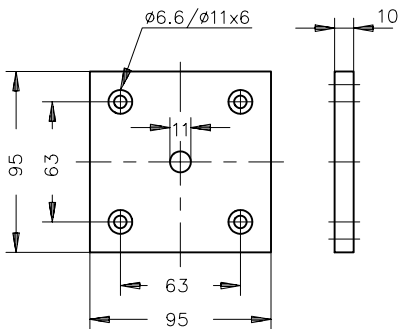
X 95 End Plate

- Caps ends of profiles / acts as a stop for carriers
- For mounting X 95 profile structures on legs or vibration isolators
- Includes 6 ea. M6 x 16 Allen-head screws

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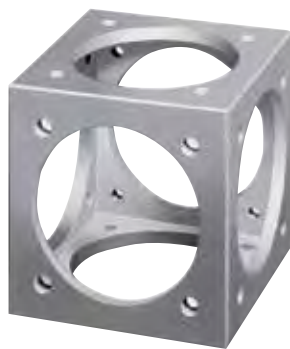X 95 Corner Connector

- Cubic coupler for assembling framework structures
- 5 faces each have diameter 6.6 mm holes for bolting coupler onto end of X 95 profile
- Other face has 4 M6 tapped holes for installing a X 95 Clamping carrier
- 1 kN load capacity
- Includes 12 ea. M6x16 mm and 4 ea. M6x25 mm screws

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X 95 Cover Plate

- For blanking off unused faces of corner connectors or open ends of X 95 profiles
- Supplied with 4 fastening screws
- Black anodized aluminum
- For volume orders have the cover plate engraved with your logo!

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X 95 Single Leg

- Single leg for X 95 rack constructions
- Height adjustable (approx. ± 10 mm)
- Requires carrier X 95 - 30 or carrier X 95 - 50 for mounting (not included, please see section Carriers)
- Swivel levelling pads supplied with screws

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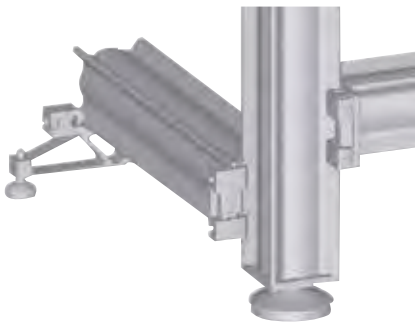
CAD files of all mechanical components can be found on the respective product sheets at www.qioptiq-shop.com below the button "Docs and Drawings".

Microbench
Nanobench
Tilt System C
Positioners

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
X 95 Double Leg

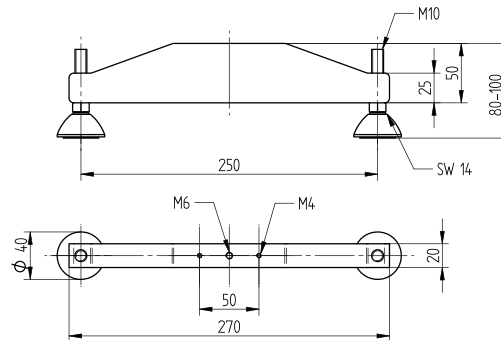- Double leg for X 95 rack constructions
- Swivel levelling pads
- Height adjustable (approx. ± 10 mm)
- Requires carrier X 95 - 30, G02 6426 000 for mounting (not included, please see section Carriers)
- Supplied with screws

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
X 95 Hinged Connector

- Allows for angle variation between two X 95 profile rails
- Used as a cross brace for X 95 structures
- $\pm 90^\circ$ pivot
- Supplied complete with mounting screws

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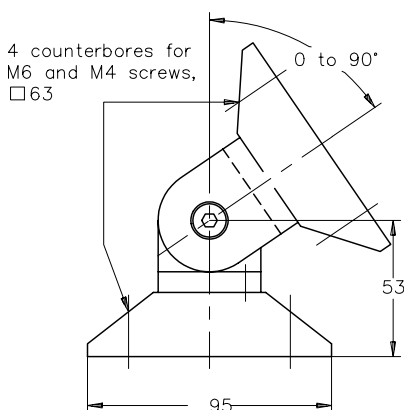Our quality criteria:

- *Stable design*
- *High-quality surface finishing*

Mounts & Posts

Qioptiq has everything you need for your optical assemblies. Our comprehensive range of mounts, posts and holders, all mutually compatible, includes an exceptionally broad range of object holders for rectangular or cylindrical construction elements, as well as optic holders for mounted and unmounted optics or prisms.

To mount optic holders on optical tables, we recommend our support rods and columns. Thanks to our large selection of base plates, clamps and magnetic feet, you can implement just about any mounting solution you want. From the clamping fork with the Easy Lift system to support rods to compatible connection clamps, Qioptiq offers a complete assortment for working with optical tables.

In addition, columns 25 and 38 are extremely sturdy mounting accessories that are ideal for heavy construction elements.

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Special Features:

A fitting gives you the option of lifting and moving the clamping fork hanging on the mount. This makes it possible to mount the clamping fork even in narrow positions in your structure or on optical tables.

Ideal Areas:

Structures on optical tables; flexible positioning and height adjustment.

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Content

Mounts and Holders compatible to FLS 95	
<i>Adjustable Holder 35-M</i>	194
<i>Adjustable Holder 55-M</i>	194
<i>XY Adjustment 50</i>	195
<i>Optic Holder 63-M</i>	195
<i>Mounting Rings for Mounted Optics</i>	196
<i>Mounting Ring for Microbench</i>	196
<i>Iris Diaphragm 50 in Mounting Ring</i>	197
<i>Plate Holder</i>	197
Mounts and Holders different heights	
<i>Centering Lens Holder 100</i>	198
<i>Object Holder 50-M</i>	198
<i>Frame Holder 100-M</i>	199
Posts, Columns and Rods	
<i>Clamping Fork</i>	199
<i>Mounting Column 14</i>	200
<i>Adjustable Height Column 75</i>	200
<i>Rods Ø 14</i>	201
<i>Connecting Clamp, 14 mm Diameter</i>	201
<i>Connection Clamp 14 V</i>	202
<i>Stop Ring, 14 mm ID</i>	202
<i>Mounts for Post Use</i>	202
<i>Post 25-200 with Clamp</i>	203
<i>Angle Bracket 40</i>	204
<i>Post 38</i>	204
<i>Stop Ring 38</i>	205
<i>Angle Bracket 80-M</i>	205
<i>Mounting Base</i>	206
Base Plates and Magnetic Base	
<i>Mounting Base 80x15</i>	207
<i>Base Plate 80x40-M</i>	207
<i>Magnetic Base</i>	207
<i>Magnetic Clamp 50-M6 (4 ea.)</i>	208
Mounting Plates	
<i>Lightweight Breadboards CleanTop® 75 Series</i>	208
<i>Mounting Plates</i>	209
<i>Mounting plates 25</i>	209
<i>Mounting Plate 220x40 mm</i>	210
<i>Mounting Plate 220x140-M</i>	210
<i>Mounting Plate 80x80</i>	210
<i>Mounting Plate 80x115-M</i>	211
<i>Mounting Plate MP 80.U</i>	211
Fasteners and Tools	
<i>Screw Set M4x8</i>	212
<i>Screw Set M4x12</i>	212
<i>Set of Screws M6x12</i>	212
<i>Set of Screws M6x16</i>	212
<i>Clamps SP 4</i>	212
<i>Set of Setscrews M6x16</i>	213
<i>Allen Wrench Set</i>	213

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
Mounts for Microbench and Nanobench

<i>Lens mounts for diameter 3 mm - 31.5 mm</i>	219
--	-----

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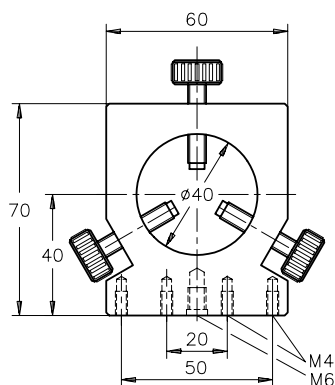
Adjustable Holder 35-M

- Holds \varnothing 10 to 35 mm radially symmetric items
- Also holds cylindrical laser heads (2 holders required)

- Nylon tipped clamping screws
- 40 mm optical axis height
- Optional grooved tips for clamping screws allow for accommodating thin-edged optics
- Mounting options: Central M6 tapped hole, M4 / 20-mm and 50-mm centers

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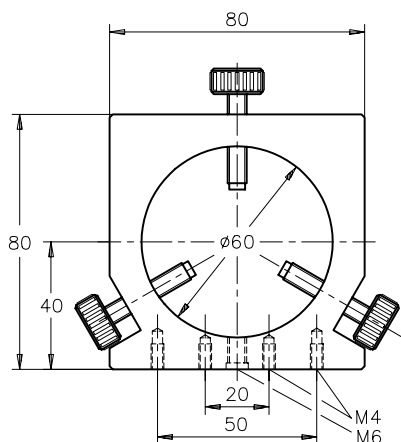Adjustable Holder 55-M

- Holds \varnothing 25 to 55 mm radially symmetric items
- Also holds cylindrical laser heads (2 holders required)

- Nylon tipped clamping screws
- Optional grooved tips for clamping screws allow for accommodating thin-edged optics
- Mounting options: Central M6 tapped hole, M4 / 20-mm and 50-mm centers

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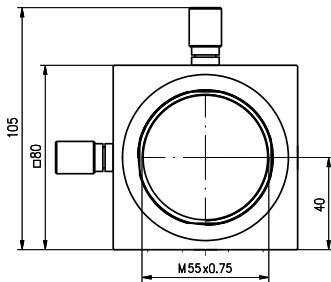
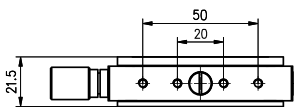
Please view chapter "Lasers and Light Sources" for appropriate lasers.

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XY Adjustment 50

The XY Adjustment 50 enables precise XY-adjustment of a wide range of optical components. For example, unmounted optics with diameters of 40 mm, 50 mm and 2" with threaded rings can be fastened directly in the adjustment. Mounts and iris diaphragms with M55 connection threads can also be adapted.

The XY Adjustment 50 can be mounted onto the rail system 65 and 95 with M4 connection threads.


The adjustment is equipped with two fine-adjustment screws with 0.25 mm pitch.

The spindle has a hexagon socket and can be driven by the provided knurled hexagon knob. If the hexagon socket is not needed, the knob can be glued in.

- Travel +/- 2.5 mm
- Sensitivity 10 µm
- Inner thread M55 x 0.75

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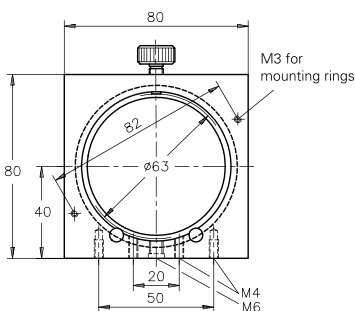
Optic Holder 63-M

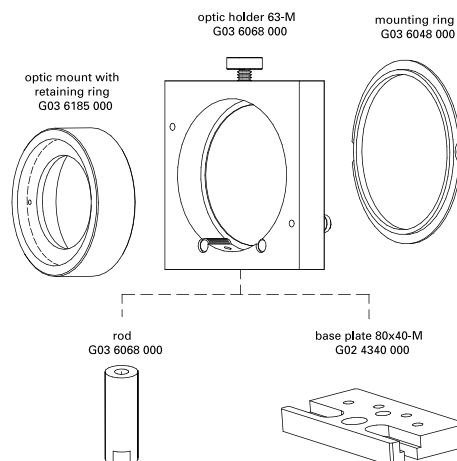
- Allows rapid interchange of optics in 63-mm outer diameter mounts
- 3-point mounting on nylon rods
- Ø 63 mm mating aperture
- 40-mm optical axis height
- Accepts unmounted optics having 5 mm or greater edge widths
- Mounting options: Central M6 tapped hole, M4 / 20-mm and 50-mm centers

Example of combinations and mounting options of optic holder 63-M

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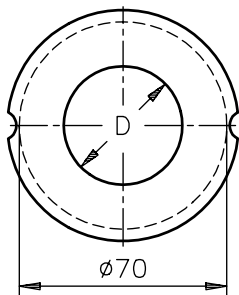
Mounting Rings for Mounted Optics

- Accepts optic mounts equipped with threaded lock rings
- The threaded holes correspond to the standardized threads of the photo industry
- To mount in Optic holder 63-M (G036068000)

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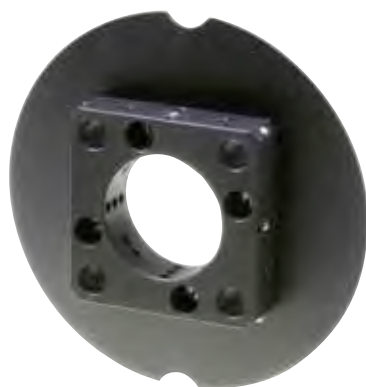
Mounting Ring for Microbench

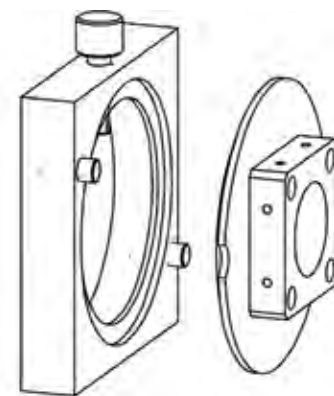
- Includes Microbench mounting plate for optic mounts with 25 mm outer diameter
- Adapts to Microbench with Microbench rods
- To mount in Optic holder 63-M (G036068000)

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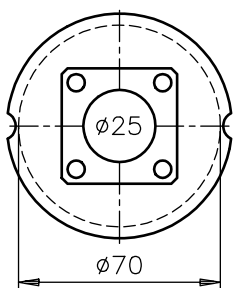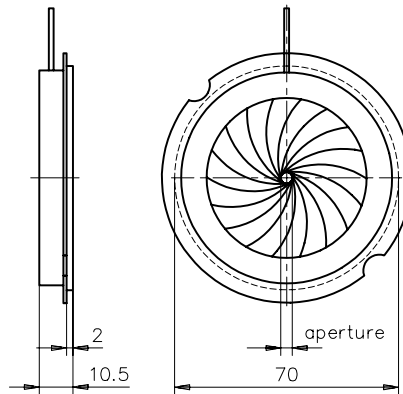
Iris Diaphragm 50 in Mounting Ring

- Opens to 50 mm diameter
- Closes down to 2.5 mm diameter
- Mates to optic holder 63-M (G036068000) and centering ring R (G070014000)

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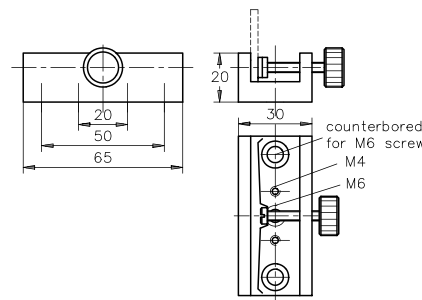
Plate Holder

- Accepts plates up to 15 mm thick
- Leaf spring for gentle clamping action
- Plastic-tipped (PA) clamping screw for rigid clamping action
- Mounting options: M6 tapped hole, counter-bored holes on 50-mm centers accept M6 screws for mounting on surfaces having M6 tapped-hole patterns, M4 tapped holes on 20-mm centers mate to Microbench mounts

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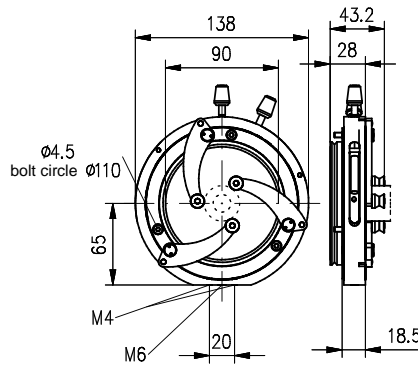
Centering Lens Holder 100

- Holds Ø 10 to 100 mm lenses
- Ø 90 mm clear aperture
- V-groove fits in adjusting ring, G07 0003 000
- Mounting options: M4 / M6 tapped hole for mounting on rods, mounting plates and carriers

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
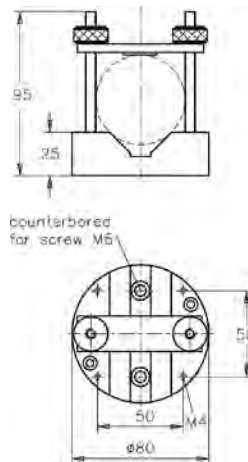
Object Holder 50-M

- Holds circular / cylindrical components having Ø 25 to 52 mm
- Accurate holder for HeNe Lasers
- Mounting options: Central M6 tapped hole, M4 / 20-mm centers carriers, Ø 6.6 / 50-mm centers

Objekthalter 50-M Object holder 50-M

Produktbezeichnung	Part No.
Item Title	Part No.
Object holder 50-M	G024446000

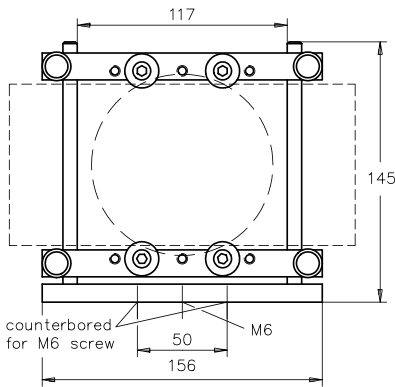Microbench
NanoBench
Tube System C
Positioners
Rail Systems

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Frame Holder 100-M

- Universal holder accepting up to 125 mm circular components or rectangular elements up to 100 mm height
- Four V-grooved rollers accommodate items having up to 8-mm edge widths
- Repositionable mounting rollers
- Base plate 30 mm wide
- Mounting options: M6 tapped hole, counter-bored holes on 50-mm centers accept M6 for mounting plates and carriers

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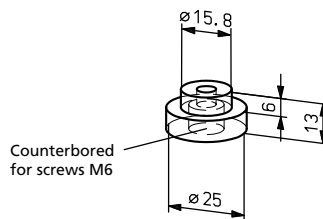
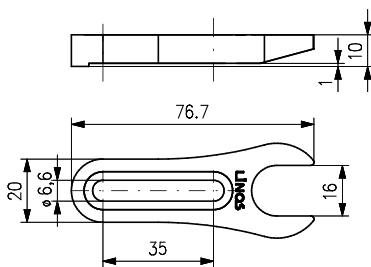
Clamping Fork

With Easy Lift System: the clamping fork can be lifted with the post, this allows an easy mounting at narrow positions on optical tables.

- For flexible mounting of components on optical tables, in connection with holder A14
- Low weight and stable
- High-strength aluminum alloy
- Clamping fork with slotted hole for M6 or 1/4" screws for attachment on optical tables
- Holder A14 suitable for M6 or 1/4" screws for attachment at posts, columns and components

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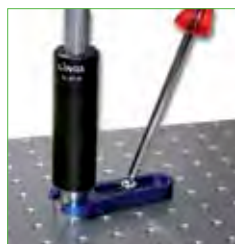
The easy lift system secures the Clamping fork on the post to avoid falling off.



Mounting of Clamping fork (with Mounting column 14 on Holder A14) on Optical Table

Easy Lift System: In case of moving the Mounting column the clamping fork clings to the column

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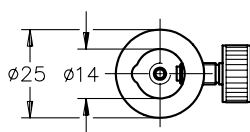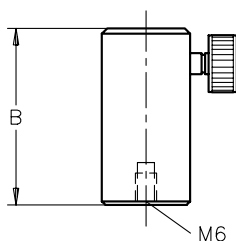
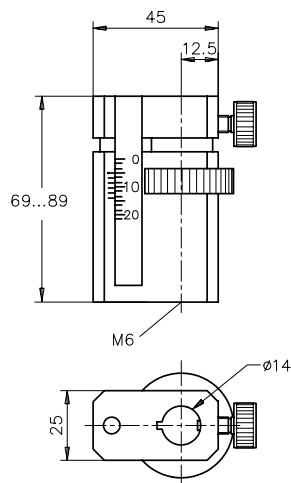
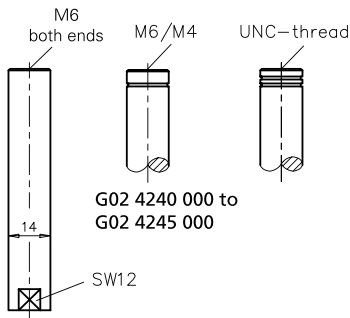
Stativstifte Ø 14 Rods Ø 14

Gewinde	Anzahl (ea.)	Länge (mm)	Part No.
Thread	Qty. (ea.)	Length (mm)	Part No.
M6/M6	1	40	G024230000
M6/M6	1	80	G024232000
M6/M6	1	120	G024234000
M6/M6	1	200	G024236000
M6/M6	1	300	G024238000
M6/M4	1	40	G024240000
M6/M4	1	80	G024242000
M6/M4	1	120	G024244000
M6/M6	10	40	G024231000
M6/M6	10	80	G024233000
M6/M6	10	120	G024235000
M6/M4	10	40	G024241000
M6/M4	10	80	G024243000

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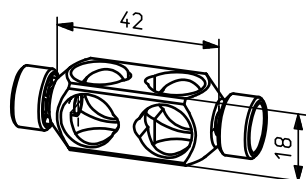Connecting Clamp, 14 mm Diameter

- For clamping Ø 14 mm rods
- Clamps rods in parallel or orthogonal orientations
- Rigid, biaxial clamping action

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
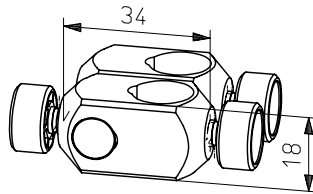
Connection Clamp 14 V

- For support rods 14 mm
- Adjustable at any angle
- Latchable adjusted angle

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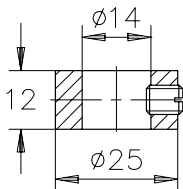Stop Ring, 14 mm ID

The stop ring makes the mounting and positioning of clamps at the rod 14 easier. On the one hand, the stop ring acts as a positioning aid, but also prevents that from inadvertently sliding components after loosening the clamping screw up on the table and get damaged.

- Adjustable stop, fits on rods with Ø 14 mm

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
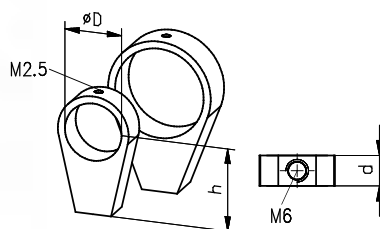
Mount 16 and Mount 25

Mounts for Post Use

These mounts allow the assembly of optics onto posts.

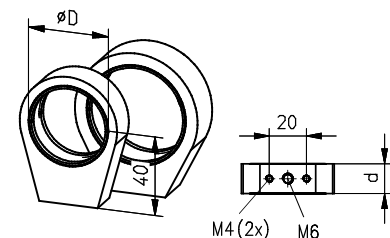- Diameter 16 mm to 35 mm for use with mounted optics of nano- or microbench mounts
- Diameter 40 mm to 80 mm for unmounted optics is delivered with retaining rings

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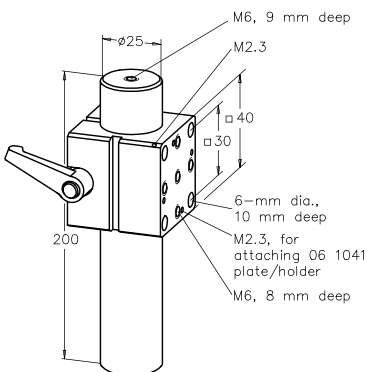
Post 25-200 with Clamp

- Mounts on optical tables or breadboards
- 25 mm diameter post, 200 mm long
- With M6 tapped hole on each end face
- Microbench rods attach directly to clamp
- M6 tapped hole pattern
- Tapped holes for attaching G061041000 mounting plate
- Cube 25 and mounting plate 25 (G061010000) attach directly to post

- Post surface is case hardened and hard-chromed
- Post surface matt, diffusely reflecting
- Mounting screws included

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
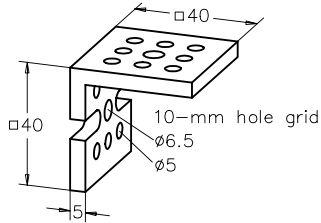
Angle Bracket 40

- Screws onto clamping fixture of post 25-200 with clamp, G061200000
- Universal bracket for assembling Microbench-compatible components

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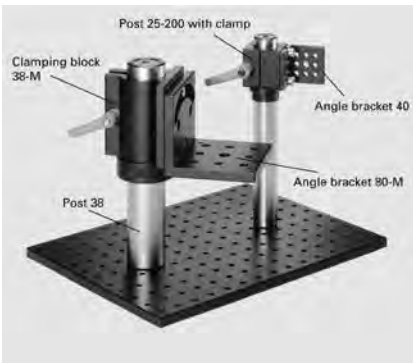
Post 38

- Solid Ø 38.1 mm steel post
- For use in combination with clamping block 38
- Metric tapped holes on both ends
- Surface case hardened to 62 HRC
- Hard-chromed matt surface
- Mounting screws included


Säule 38 Post 38

Produktbezeichnung	Gewinde	Gewicht (kg)	Höhe (mm)	Part No.
Item Title	Thread	Weight (kg)	Height (mm)	Part No.
Post 38-200-M	M6	1.76	200	G024944000
Post 38-350-M	M6	3.1	350	G024946000

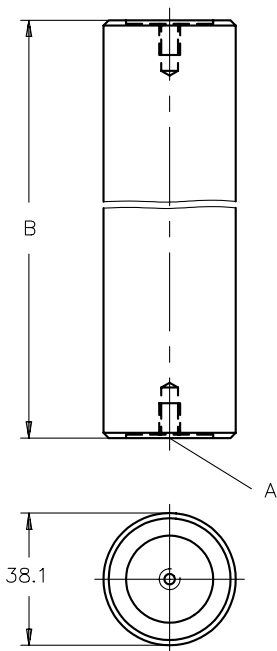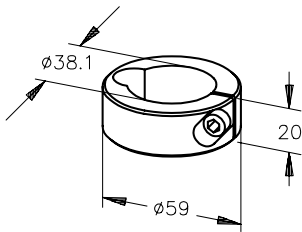Microbench
nanobench
Tube System C
Positioners
Rail Systems

Stelling 38

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Stop Ring 38

- Firmly clamps on post 38 and acts as stop

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

Angle Bracket 80-M

- Universal angle bracket
- Mates to optical tables, and flat rail carriers
- Components mount from either side

Montagewinkel 80-M Angle bracket 80-M

Produktbezeichnung	Part No.
Item Title	Part No.
Angle bracket 80-M	G024334000

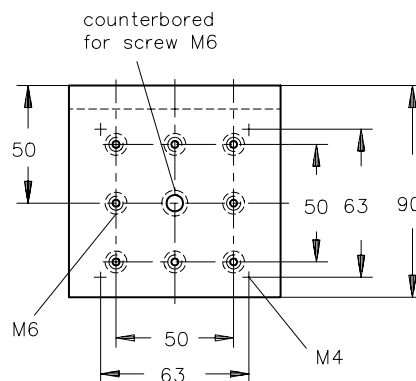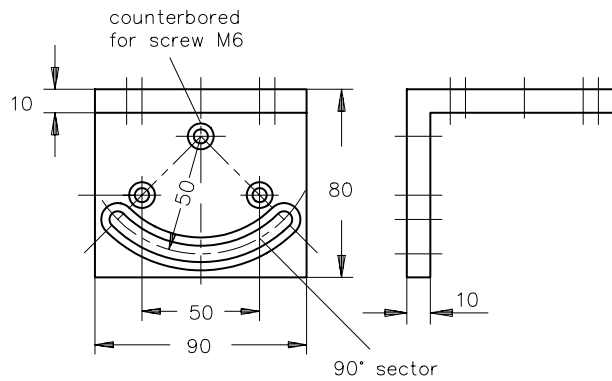Application example for mounting base

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

Mounting Base

- Stable base for mounting posts (such as post 38) or vertically mounting X 95 profiles
- Machined flat mounting surface, 95 mm wide
- M6 threaded mounting holes for mounting plates and low profile rail, G022080000 etc.
- Levelling feet with a height adjustment up to 20 mm

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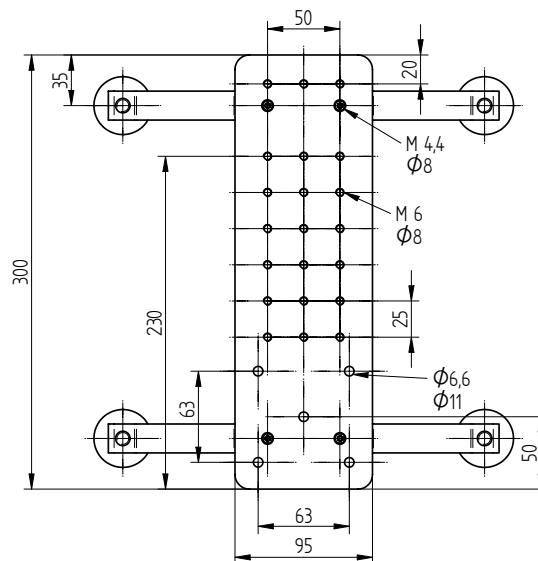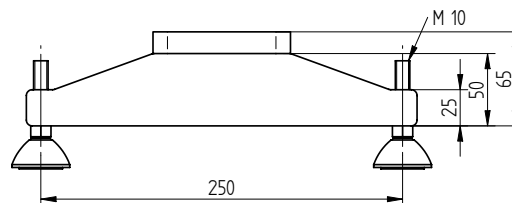
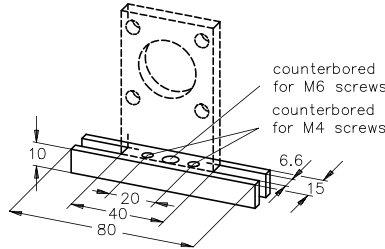
Mounting Base 80x15

- For mounting Microbench system components on optical tables
- Allows narrower spacings between components

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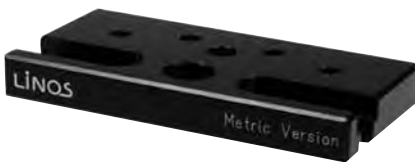
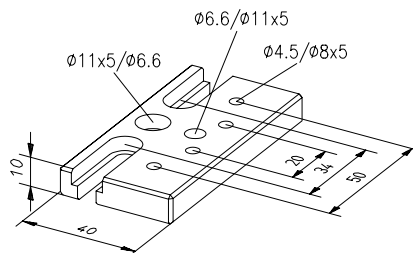
Base Plate 80x40-M

- Universal base plate accepting narrow components
- Tapped holes accept Microbench components and components with 50 mm mounting grid
- Mounting screws included

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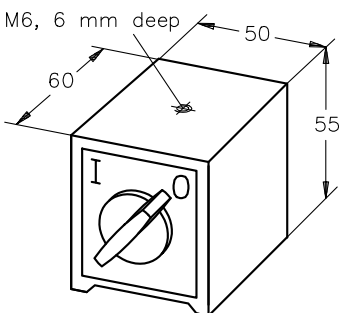Magnetic Base

The Magnetic Base can be switched on and off as needed. With the magnetic current activated, any of a broad range of assembly components can be firmly attached to a ferromagnetic surface. When the magnetic base is switched off, the magnetic current is cut and the base can be moved effortlessly into another position. When the magnetic current is on, it has a holding force of 800 N to ensure high stability. The base has a centered M6 tapped bore to attach columns for support rods or other components.

- Accepts columns, mounting plates, etc.
- Equipped with magnet-release knob
- Approx. 800 N gripping force
- Central M6 tapped hole

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
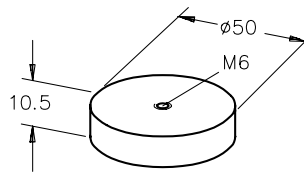
Magnetic Clamp 50-M6 (4 ea.)

- For posts, mounting plates etc.
- Multiple use produces high gripping force
- Rubber cuff (exchangeable) for higher shear force and protection of working surface
- Gripping force of 200 N per magnet (without rubber cuff)
- Mounting screws and cupped washers included

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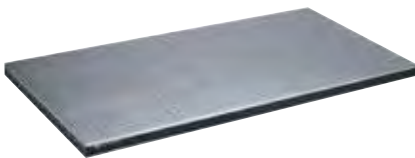
Lightweight Breadboards CleanTop® 75 Series

Minimizing the weight of Optical Tables is becoming more of a consideration for many applications. The development of new manufacturing techniques has resulted in a new, ultralight design.

- TMC breadboards in ultra-light design
- Steel honeycomb core
- Top and bottom plates made of stainless steel
- Countersunk tapped bores with 3 thread turns
- Economically priced
- Core: Steel honeycomb, closed cell, 0.2 mm thick foil
- Core shear modulus: 19,300 kg / cm²
- Core density: 230 kg / m³
- Flatness: ±0.13 mm
- Top skin: 430 series ferromagnetic stainless steel, 2 mm thick
- Bottom skin: 400 series ferromagnetic stainless steel, 2 mm thick
- Sidewalls: High-pressure laminate
- Tapped holes: M6, backed by CleanTop nylon cups, 3 threads

Leicht-Breadboards CleanTop® Serie 75 Lightweight Breadboards CleanTop® 75 Series

Produktbezeichnung	Gewicht (kg)	Part No.
Item Title	Weight (kg)	Part No.
Leicht-Breadboard 450x600x50mm, M6	14	G437510312
Leicht-Breadboard 450x1200x50mm, M6	28	G437510412
Leicht-Breadboard 600x600x50mm, M6	17	G437511312
Leicht-Breadboard 600x900x50 mm, M6	25	G437511512
Leicht-Breadboard 600x1200x50mm, M6	34	G437511912
Leicht-Breadboard 900x1200x50mm, M6	50	G437513512

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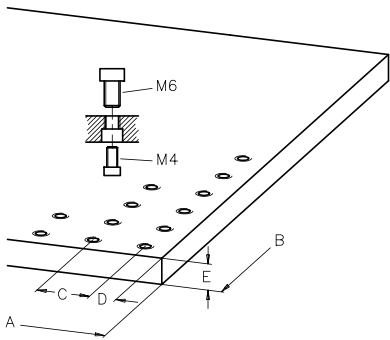
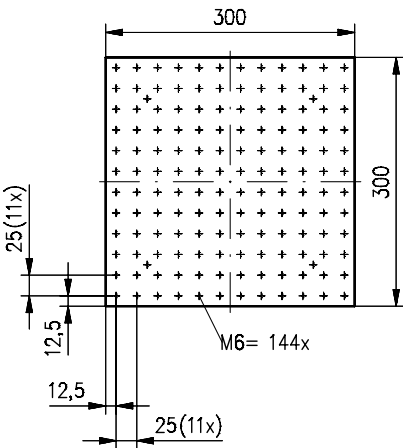
Mounting plates 25


- *Stable mounting plates for compact setups*
- *M6 x 25 tapped hole pattern*
- *Thickness 10 mm*
- *Non-slip rubber feet included*
- *Solid aluminum construction*
- *Black anodized*

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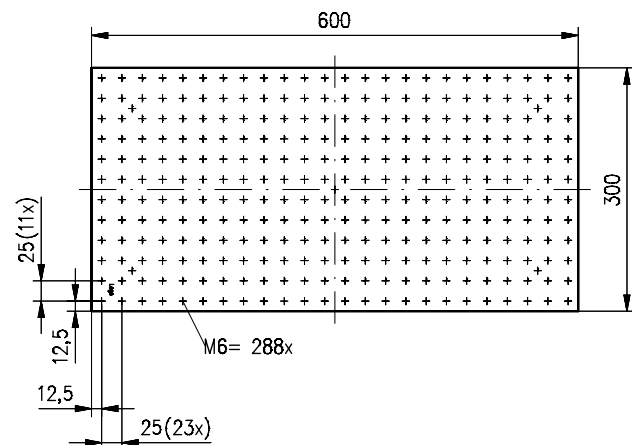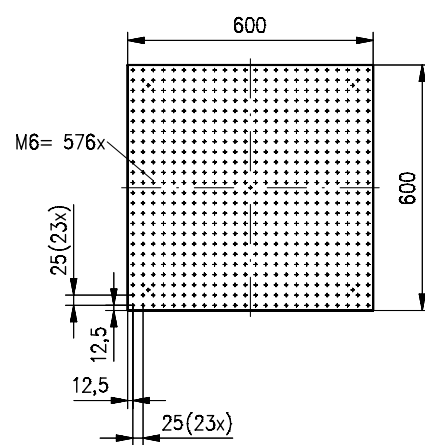
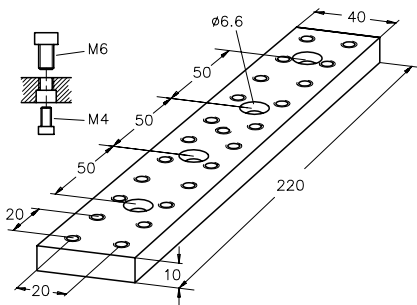
Mounting Plate 220x40 mm

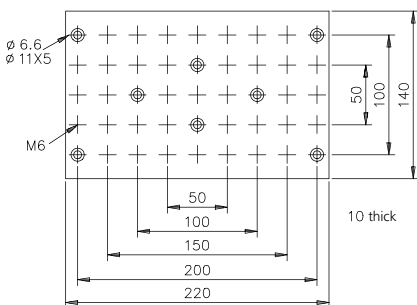- Used as a base plate for Microbench structures
- Mounts directly to carriers
- 20 mm hole pattern
- Top threaded M6, bottom counter bored for M4
- Used as a base plate for linear translation stages G065114000 and G065115000

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
Mounting Plate 220x140-M

- For assembling sets of components and modular structures
- Fit to post 38
- May be used as separate table surface

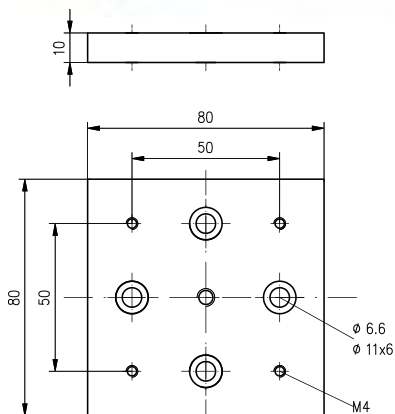- Grid 25 mm
- See section Posts, Columns and Rods regarding post 38

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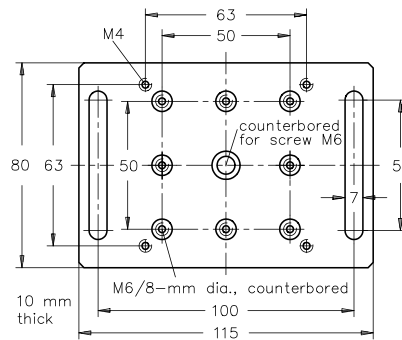
Mounting Plate 80x80

- Mounting plate to attach manual positioners with hole pattern 50 mm M4 to optical tables or X95 carriers with M6 at pattern 25 mm

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Mounting Plate 80x115-M

- Universal mounting plates equipped with slots
- Metric / inch tapped-hole patterns
- Components mount from either side using tapped holes provided / allen-head screws

Montageplatte 80x115M Mounting plate 80x115-M

Produktbezeichnung	Part No.
Item Title	Part No.
Mounting plate 80x115-M	G024330000

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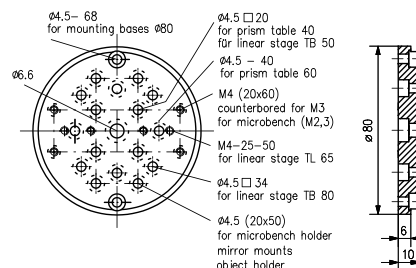
- Ø 80 mm
- Dicke 10 mm

Mounting Plate MP 80.U

- Adapter plate suitable for diverse Microbench components e.g. holder 25 / 30 / 35, rod holder, bench plates
- For prism table 40S/60S
- For linear stages TB 32 / TB 50 / TB 80
- For diverse object and mirror holders
- Mountable on mounting bases without clamps
- To assemble on carrier FLR 95 and carrier X 95 with help of clamps SP 4, G085922000


- Ø 80 mm
- 10 mm thick

Montageplatte MP 80.U Mounting plate MP 80.U

Produktbezeichnung	Part No.
Item Title	Part No.
Mounting plate MP 80.U	G024988000

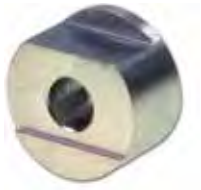Example: Mounting plate on X95 carrier

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
Clamps SP 4

The clamps can be used to mount round mounting plates with a diameter of 80 mm onto carriers having suitable tapped-hole patterns M4 square 63 mm. The mounting plates can be rotated to the desired position.

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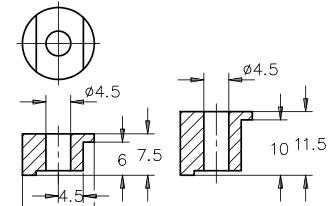Screw Set M4x8

- Consisting of:
- Hex socket head screws M4x8 (ISO7380), 50 ea.
- Cupped washers D8/D4.0 (DIN433), 50 ea.
- Wrench 2.5 mm, 1 ea.

Produktbezeichnung	Part No.
Item Title	Part No.
Screw set M4x8	G061255000

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Screw Set M4x12

- Consisting of:
- Hex socket head screws M4x12 (ISO7380), 100 ea.
- Cupped washers D8/D4.0 (DIN433), 100 ea.
- Wrench 2.5 mm, 1 ea.

Produktbezeichnung	Part No.
Item Title	Part No.
Screw set M4x12	G061256000

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Set of Screws M6x12

- Consisting of:
- Hex socket head screws M6x12 (ISO7380), 100 ea.
- Cupped washers D11/D6,3 (DIN433), 100 ea.
- Wrench 4 mm, 1 ea.

Produktbezeichnung	Part No.
Item Title	Part No.
Set of screws M6x12	G026665000

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Set of Screws M6x16

- Consisting of:
- Hex socket head screws M6x16 (ISO7380), 50 ea.
- Cupped washers D11/D6,3 (DIN433), 50 ea.
- Wrench 4 mm, 1 ea.

Produktbezeichnung	Part No.
Item Title	Part No.
Set of screws M6x16	G026666000

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Set of Setscrews M6x16

- Consisting of:
- Setscrews M6x16 (DIN913), 50 ea.
- Wrench 3 mm, 1 ea.

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Allen Wrench Set

- Consists of:
- Wrenches of the sizes 1.5; 2; 2.5; 3; 4; 5; 6; 8; 10

- Chrome-vanadium, nickel-plated
- In plastic clip with hanger

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Früher
In former times

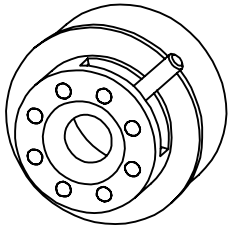
Montage und Mechanikfertigung in
 Göttingen um 1910
*Mounting and production of mechanical
 components in Goettingen about 1910*

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Iris Diaphragm Nanobench

- Min. aperture 0.5 mm, max. aperture 7 mm
- Outer-Ø 16 mm to insert in Mounting plate N 16, Holder N 40 and Cube N

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

Iris Diaphragms Microbench


- Adjusting lever screwed-in (M1.6)
- Outer diameter 25 mm
- G061657000 with completely closing aperture allows use as dimmer and switch

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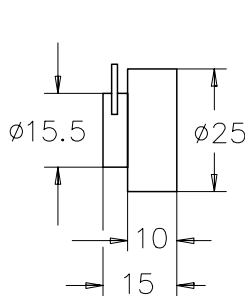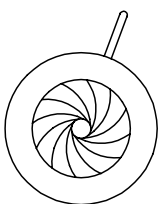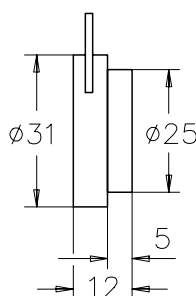
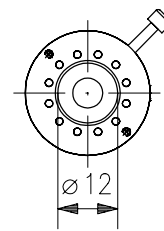
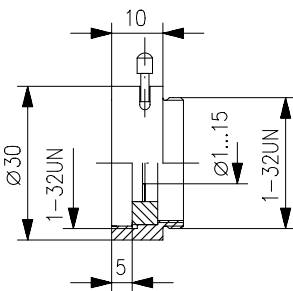
Iris Diaphragm Tube Mounting System C


- Free aperture \varnothing 1 - 15 mm or \varnothing 1.2 - 22 mm
- Adjusting lever screwed-in (M1.6)
- Outer \varnothing 30 mm or 35 mm
- Internal and external 1-32UN (C Mount) or internal M32.5 x 0.75 on both sides (G061653000)

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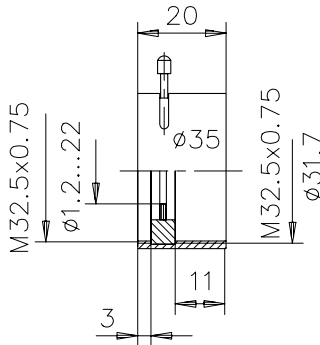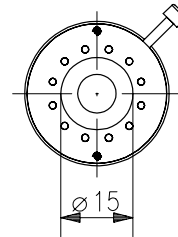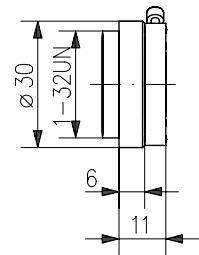
*) Completely closing iris diaphragm: It is practical to use completely closing iris diaphragms on an optical bench, for instance, if you need to completely block off a laser beam path. Iris diaphragms are also ideal for use as a combined dimmer and switch for lasers or advanced arc lamps, which cannot be briefly switched on or off once they are triggered.

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

Iris Diaphragms, Completely Closing

It is practical to use completely closing iris diaphragms if you need to completely block off a laser beam path. Iris diaphragms are also ideal for use as a combined dimmer and switch for lasers or advanced arc lamps, which cannot be briefly switched on or off once they are triggered.

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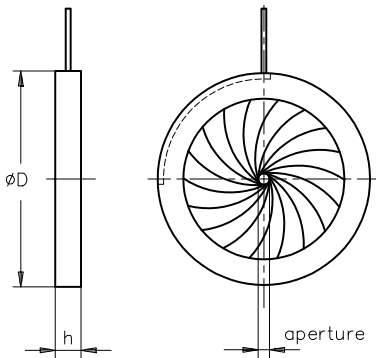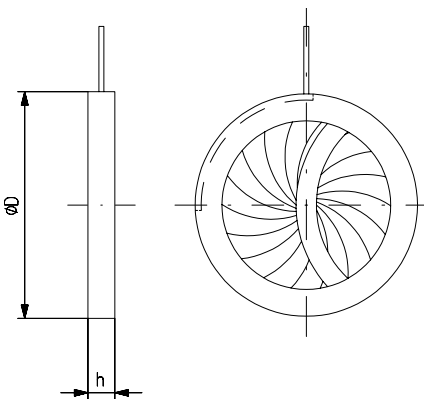
Iris Diaphragms with Different Outer Diameters

- Outer diameter from 24 mm up to 120 mm
- Min. aperture from 1.0 mm up to 3.5 mm
- Including adjusting lever

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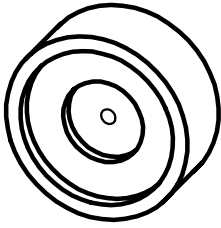Mounted Pinholes

- For diffraction experiments, alignment purposes, projection applications
- Supplied in \varnothing 25 mm mounts for Microbench and \varnothing 16 mm mounts for Nanobench
- Chemically etched apertures
- Chemically blackened brass
- 0.1 mm plate thickness
- Coaxiality tolerance 0.1 mm

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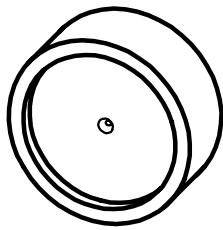
For adjustment in Microbench or Nanobench system the Centering Mounting Plate (G06 1025 000, chapter Microbench) and XY Translation stage N (G05 0564 000, chapter Nanobench) are ideally suited.

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
Pinholes for Spatial Filters

- For spatial filters
- For spatially filtering laser beams
- Reflex free types available

- Supplied in \varnothing 25 mm mounts for Microbench or in \varnothing 16 mm mounts for Nanobench
- Fabricated from platinum-iridium or molybdenum

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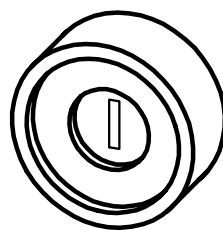
Slits

- For diffraction experiments, alignment purposes, projection applications

- Supplied in \varnothing 25 mm mounts for Microbench or in \varnothing 16 mm mounts for Nanobench
- Chemically etched slits
- Chemically blackened brass
- 0.1 mm plate thickness
- 5 mm height of slit

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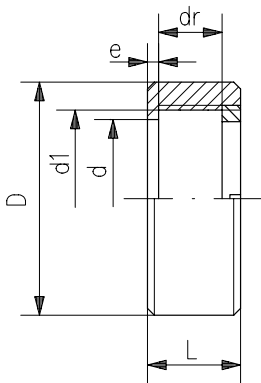
Lens mounts for diameter 3 mm - 31.5 mm

These lens mounts are ideal for mounting optics – whether from LINOS or from other manufacturers – with diameters of 3 mm to 31.5 mm. A threaded ring for fastening the optics is included with the mount. These lens mounts make it easy to integrate optics in LINOS Microbench and Nanobench systems, as well as in holders for mounting on support rods.

- Suitable for the Nanobench (N) and Microbench (CL for lenses, CA for achromats) system
- Retaining ring included
- Tip: LINOS optics can be mounted before shipping. In this case, the optics designation is engraved on the mount, which can prevent confusion. Pre-mounting also protects the optics from the damage that can occur when switching lenses in a mount.

Optikfassungen Durchmesser 3 mm - 31.5 mm Lens mounts for diameter 3 mm - 31.5 mm

Produktbezeichnung	d (mm)	E (mm)	Ø D (mm)	Ø d1 (mm)	L (mm)	dr min. (mm)	dr max. (mm)	Part No.
Item Title	d (mm)	E (mm)	Ø D (mm)	Ø d ₁ (mm)	L (mm)	dr min. (mm)	dr max. (mm)	Part No.
Mount N3	2.5	0.5	16	3	8	0.5	3.4	G050423000
Mount N5	4	0.5	16	5	8	0.8	3.4	G050425000
Mount N6	5	0.5	16	6	8	0.8	3.4	G050426000
Mount N8	7	0.5	16	8	8	0.8	3.4	G050428000
Mount N10	9	0.5	16	10	8	1.1	3.4	G050430000
Mount N12.5	11.5	0.5	16	12.5	8	1.1	5.9	G050432000
Mount N12.7	11.5	0.5	16	12.7	8	1.1	5.9	G050433000
Mount CL 18	17	1	25	18	10	1	4	G063653000
Mount CL 19	17	1	25	19	10	1	4	G063661000
Mount CL 22.4	21.4	1	25	22.4	10	1	6	G063651000
Mount CL 22.4	21.4	1	25	22.4	18.5	1	15	G063671000
Mount CA 18	17	1	25	18	10	3	6	G063658000
Mount CL 25.4	24	1.5	30	25.4	12	1.5	9	G063663000
Mount CL 31.5	30	3.5	35	31.5	16	1	10	G063668000

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
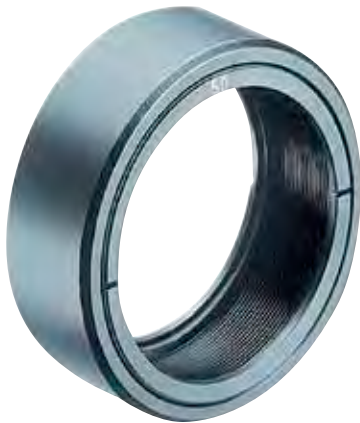
Optic Mount 63 (40, 50, 50.8 mm)

- Choice of three models for mounting Ø 40; 50; 50.8 mm optics
- Full width internal thread
- Accepts optics having 1.5 to 20 mm edge widths
- Fits optic holder 63-M, G036068000
- Supplied complete with two retaining rings

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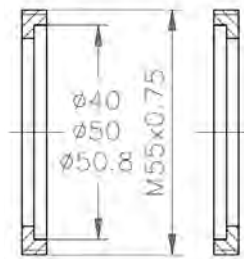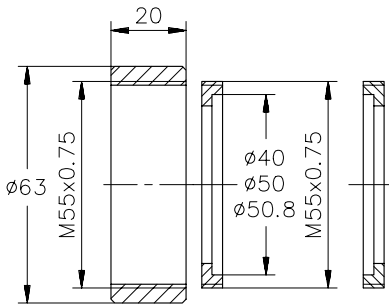
Microbench
Handbench
Tube System C
Positioners
Rail Systems

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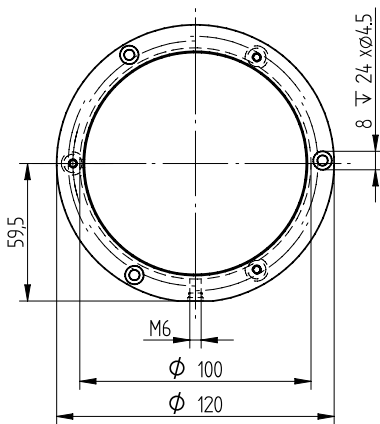Mount 100


- For mounting \varnothing 100 mm optics
- \varnothing 97 mm clear aperture
- Mounting options: Central M6 tapped hole for mounting plates or rods

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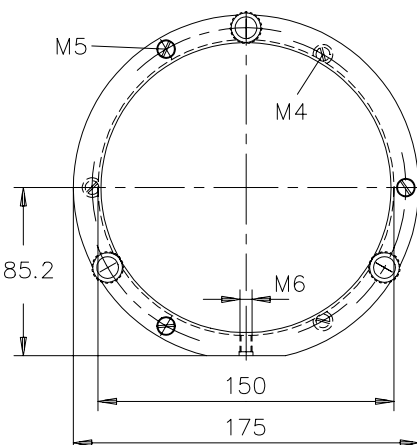Holder 150

- For mounting \varnothing 150 mm optics
- \varnothing 147 mm clear aperture
- For max. 10 mm or 28 mm optics edge widths
- Holder 150.28 inclusive adjustment screws and spring fastener
- Mounting options: Central M6 tapped hole for mounting plates or rods

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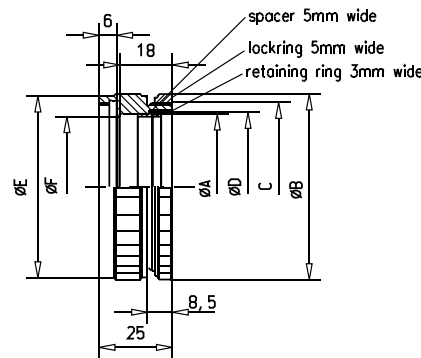
Lens Mount


- For quick and easy realization of optical systems as prototypes, models or small series
- Integration of unmounted optics with diameters 40 mm, 50 mm, 63 mm, 80 mm
- Compatible with Tube Mounting System C components by using adapters (diameter 31.5 mm, 25.4 mm, 22.4 mm and smaller)
- Possibility to connect filters, diaphragms etc. from other sources
- Connecting thread of the standard line of objectives according to DIN 13, DIN 19002 and ISO 1948 of photo industry

- Equipped with retaining ring and locking collar

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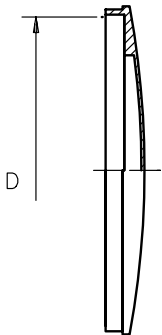Lens Protection Cap

- Protect optics against dust and damage
- Type name is equivalent to inner diameter D
- Suitable for lens mounts, adapters and iris diaphragms

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Our quality criteria:

- *High-quality adjustment screws with virtually no backlash*
- *Stable design*
- *Precise positioning*
- *Upgradeable*

LINOS Mirror Mounts

Mirror mounts are an essential element in optical assemblies. The variety of mounts available from Qioptiq meets a broad range of requirements in science and industry.

The LINOS adjust.x series is both economical and precise. Models with integrated base plates have dimensions compatible with those of Lees mirror mounts, so you can upgrade to Lees at any time. Our standard M series includes extremely compact mirror mounts handled from above, gimbal mirror mounts, and rugged industry mirror mounts with incomparable long-term stability.

Lees mirror mounts, the premium class mounts from LINOS, have a unique design that guarantees the highest possible precision. Their excellent quality is further ensured by features such as fine adjustment screws that are specially ground for mutual fit with the least possible backlash, counter surfaces of sapphire and the special Krytox lubricant.

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Special Features:

Qioptiq offers a comprehensive selection of mirrors that fit these mounts, from metallic mirrors to dielectric multi-coated mirrors.

Ideal Areas:

Position adjustment for mirrors of 0.5" to 2" in diameter; experimental setups using optical tables; installation in instruments.

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Unique aluminum base design

The solid LM1 optic mount base and plate are machined from aluminum alloy and are protected with a non-shiny black aluminum oxide coating known as anodize. The unique design of the base is more rigid and has a higher frequency than flat mounts. The LM1's wide bottom provides superior mounting to tables or accessories. Thus, greater stability with less susceptibility to vibration.

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The LM1 is equipped with three 10-80 adjustment screws each with 0.175 in. of travel minimum. The adjustment screws are considered to have the finest resolution available today.

Dynamic range

The LM1 mount can be angularly adjusted to more than $\pm 25.0^\circ$. The front plate can also translate more than ± 0.175 in.

Sapphire hard seat technology

We use a large $3/16 \times 0.034$ in. sapphire flat polished to optical standards on all LM1 mounts. The hard cylindrical vee is also standard on all LM1 mounts. These seats are so tough, they have been known to work flawlessly even when the mount has been dropped.

Nylon tipped set screws

Kinematically locates the optics and secures it without stressing or damaging it. They also allow for easy removal and replacement.

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Metric

The standard mounting hole in the base is M6. The top set screw in the front plate is 4-40 UNC or 6-32 UNC.

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LM2 Optic Mount, Compact

Compact Precision Mirror Mount for 1" optics

Unique aluminum base design

The solid LM2 optic mount base and plate are machined from aluminum alloy and are protected with a non-shiny black aluminum oxide coating known as anodize. The unique design of the base is more rigid and has a higher frequency than flat mounts. The LM2's wide bottom provides superior mounting to tables or accessories. Thus, greater stability with less susceptibility to vibration.

Stainless steel adjustment screws in phosphor bronze bushings

The LM2 is equipped with three 1/4-80 adjustment screws each with 0.5 in. of travel. The adjustment screws are considered to have the finest resolution available today.

Dynamic range

The LM2 mount can be angularly adjusted to more than $\pm 16.5^\circ$. The front plate can also translate more than ± 0.2 in.

Sapphire hard seat technology

We use a large $1/4 \times 0.031$ in. sapphire flat polished to optical standards on all LM2 mounts. The hard cylindrical vee is also standard on all LM2 mounts. These seats are so tough, they have been known to work flawlessly even when the mount has been dropped.

Nylon tipped set screws

Kinematically locates the optic and secures it without stressing or damaging it. They also allow for easy removal and replacement.

1/4-20 or M6 mounting hole

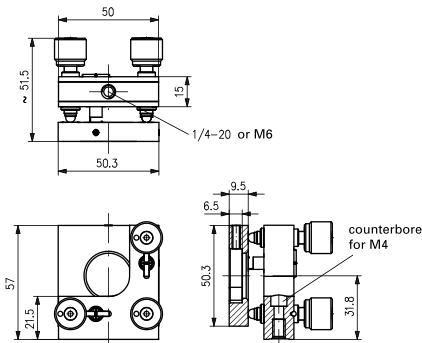
We provide the largest and strongest mounting holes to ensure stability. Deflections are solely the result of material elasticity.

Krytox fluorinated lubricant

Krytox greases have vapor pressures of 8×10^{-8} to 6×10^{-9} torr at 30°C (100°F) according to Du Pont and thus are unlikely to outgas on your expensive optics. Krytox also has a high degree of chemical inertness and outstanding oxidation resistance. Therefore, the adjustment screws will maintain their original smooth feeling if used daily, monthly, or yearly.

Metric


The standard mounting hole in the base is M6. The top set screw in the front plate is 8-32 UNC.

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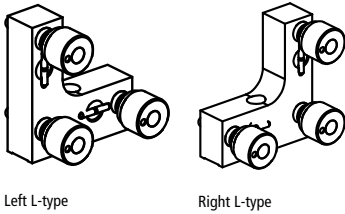Front plate, Plain without hole

Front plate for 1.0" Optic

Front plate for 1.0" Optic; Offset-Left or Offset-Right

Leas

Quick Change Indexer on Mounting Plate

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Quick Change Indexer LI2

The quick change indexer LI2 permits simple, reproducible exchange of mirror mounts or frame or plate holders in laser beams. Its catch design using three spheres ensures clean reproducibility.

The upper part can be securely offset from the lower part in 15° steps just by pressing the two pushbuttons, without the need for any extra tools. Usually, an LM2 mirror mount is attached to the upper part of the indexer so that this provides a beam height of 4".

The lower part of the indexer remains firmly attached to the optical table. Qioptiq offers quick riser bases in 2 versions: with lower base and mounting plate. Each version is available in inch or metric dimensions.

The upper and lower parts (upper and lower base parts) can also be separately supplied so that several mirror mounts can be used in exact positions on a lower part or one mirror mount can be attached in different positions. This is advantageous, if, for instance, mirrors with different coatings need to be used in succession.

The LI2 indexer can be installed in a vertical, horizontal or even in an overhead position, guaranteeing the best possible stability and reproducibility in any installation position. The mounting plate is designed to accommodate spacing of 25.4 mm or 25 mm, which is commonly used for most optical tables.

The indexer enables X-Y translation using the slots in the base plate and convenient and secure attachment by an excenter fixing. The indexer is a strictly mechanical device for mounting optical components and does not have any magnet. The LI2 is lubricated with Krytox to prevent outgassing.

- Reproducible interchange of mirror mounts
- Ideal alternative to flip mirror mounts
- Easy handling

Microbench
nanobench
Tilt System C
Positioners
Rail Systems
Mounts and Posts

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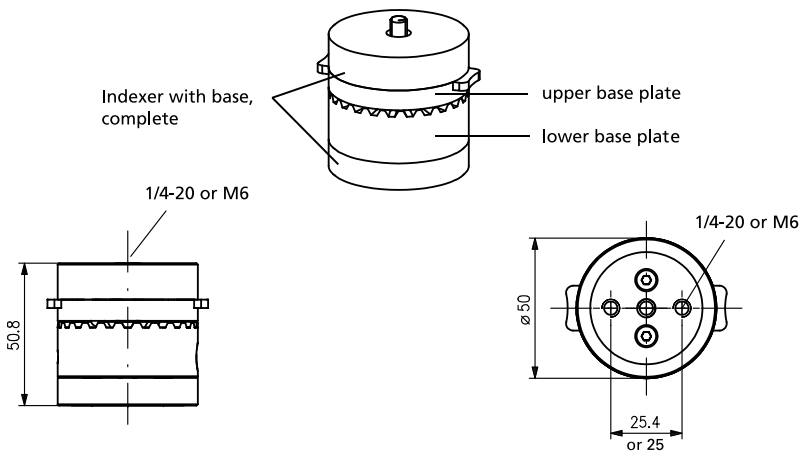
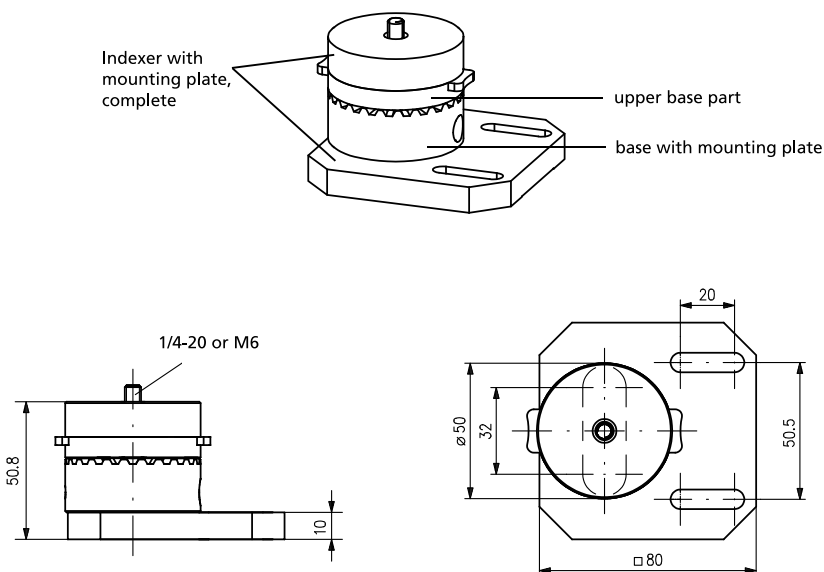
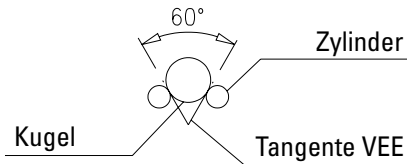
Hard Seats

Lees Optical Instrument Co., Inc. was the first to offer sapphire hard seat technology when the hard seat option was developed for LM1 and LM2 optical mounts. This technology is standard on all LM1, LM2, and LM3 optical mounts.

Sapphire technology was chosen for its hardness, second only to diamond, and an optical quality polished surface finish. The sapphire disk is available in 3/16 or 1/4 inch diameters. The vee is also unique in that it is constructed from two hardend and lapped steel cylinders. The cylinders are arranged to form a 60° vee at the tangency point of the ball and cylinders.


Sapphire technology advantages:

- Extremely hard
- Highly polished surface
- Low friction
- Smooth bidirectional translation


Hard vee advantages:

- Very hard
- Lapped surfaces
- Precision vee form
- Precision translation
- Low friction
- Incorporated in brass-base

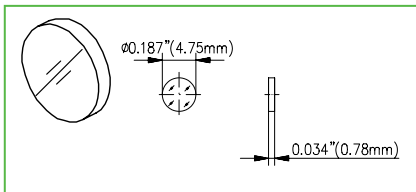
HS010-10


HS250-10

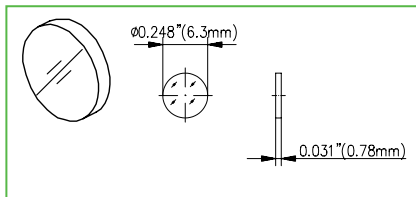
HS010-20

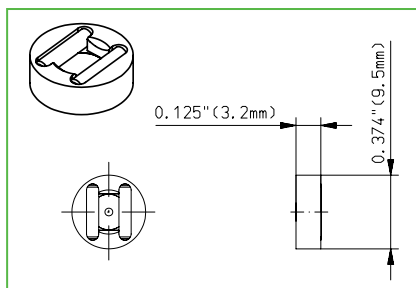Ball LM1

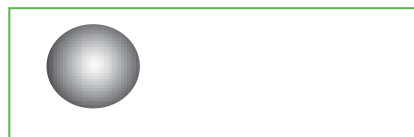Ball LM2-3

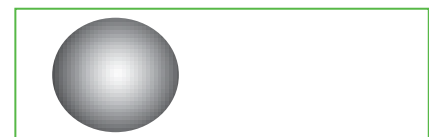Gegenlager Hard Seats


Produktbezeichnung	Part No.
Item Title	Part No.
Sapphire hard seat HS010-10	G404600000
Sapphire hard seat HS250-10	G404601000
Ball cell HS010-20	G404602000
1/8 inch ball LM1; 5 ea.	G404604000
3/16 inch ball LM2-3; 5 ea.	G404605000

Microbench
Nanobench
Tube System C
Positioners
Rail Systems
Mounts and Posts

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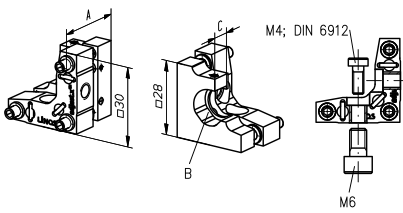adjust.x Small L-Type

Mirror mounts in the adjust.x series are available in three different designs to suit individual requirements.
 The adjust.x small mount has an L-shaped base with M6 connection threads and counterbores for M4 screws. It can be fastened over both of the brackets on the base.
 The mirror mounts are equipped with three fine adjustment screws with 0.15 mm thread pitch Allen screws.

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- 1: Closed front plate for 1/2" mirror. The mirror is locked in place by a threaded pin with nylon tip.
- 2: Closed mounting plate for 1" mirror. The mirror is locked in place by a threaded pin with nylon tip.
- 3: Open mounting plate for 1/2" mirror. The open design lets you direct a beam past a mirror at extremely close proximity. Also suitable for beam-splitter plates.

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adjust.x Medium L-Type

Mirror mounts in the adjust.x series are available in three different designs to suit individual requirements.

The adjust.x medium mount has an L-shaped base with M6 connection threads and counterbores for M4 screws.

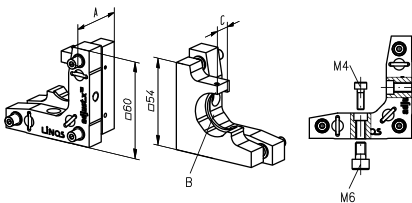
It can be fastened over both of the brackets on the base. The mirror mounts are equipped with three fine adjustment screws with 0.15 mm thread pitch allen screws.

Three types of front plate are available:

1: Closed front plate for 1" mirror. The mirror is locked in place by a threaded pin with nylon tip.

2: Closed mounting plate for 2" mirror. The mirror is locked in place by a threaded pin with nylon tip.

3: Open mounting plate for 1" mirror. The open design lets you direct a beam past a mirror at extremely close proximity. Also suitable for beam-splitter plates.

	A (mm)	B (mm)	C (mm)
adjust.x LT m 1"	36.5	25.4 (1")	10
adjust.x LT m 1" RR	36.5	25.4 (1")	3-8.5
adjust.x LT m 1" OF	36.5	25.4 (1")	10

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
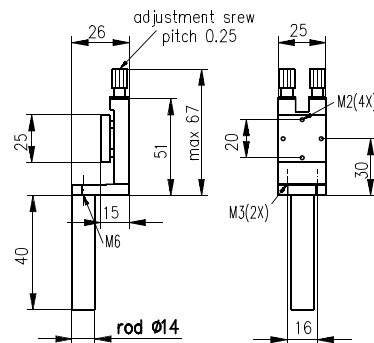
Mirror Mount MF-25PF

- Top mounted adjustment screws for space restricted setups
- Two-axis tilt control
- Kinematic Movement
- 25 x 25 mm mounting area
- Precision adjustment screws with 0.25 mm pitch
- Rod Ø 14 mm included

**Spiegelhalter MF-25PF
Mirror mount MF-25PF**

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Mirror mount MF-25PF	G404195000

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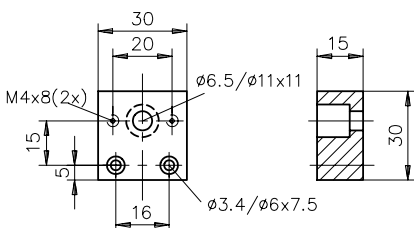Adapter Plate MF-25

- To raise the beam height to 40 mm on direct mounting the Mirror mount MD-25FP or MF-25PF on carriers and mounting plates

**Adapterplatte MF-25
Adapter plate MF-25**

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Adapter plate MF-25	G404201000

Mirror and Mounting column 14 not included

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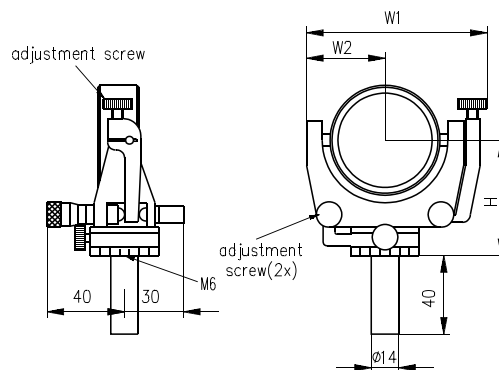
Gimbal Mirror Holder MA

- Precision micrometer movement
- Coarse angular translation, 360°
- Fine angular translation, 0.25 mm pitch
- Color coded knobs
- Rod $\varnothing 14$ mm included

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These mirror mounts offer the advantage that the surface of the mirror in the center point is always at the rotation point. This means adjustment avoids translation movement of the rotation point.

Gimbal-Spiegelhalter MA Gimbal Mirror holder MA

Produktbezeichnung	$\varnothing D$ (mm)	W1 (mm)	W2 (mm)	Max. Spiegeldicke (mm)	H (mm)	Part No.
Item Title	$\varnothing D$ (mm)	W1 (mm)	W2 (mm)	Max. mirror thickness (mm)	H (mm)	Part No.
Gimbal mirror holder MA	25.4	71	30	6	47	G404211000

Microbench
NanoBench
Tube System C
Positioners
Rail Systems
Mounts and Posts

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Our quality criteria:

- *Reliable: performance tested in the lab*
- *Compatible: can be combined with established LINOS catalog components, especially with Microbench systems*
- *Fast: complete assembly kit for standard lab applications*
- *Beneficial: inexpensive due to set-price*

LINOS Q-Sets

A Q-Set is an assembly of components from the LINOS catalog for specific lab applications. Your benefit: Save time since we select the required components from the catalog portfolio for your optical or optomechanical system for you. Easy assebmling is guaranteed!

Each Q-Set is tested thoroughly. By that we assure the high reliability of Qioptiq components also for this product line. To meet preferably all requirements and to address a broad spectrum of applicants a Q-Set is usually designed for a broad wavelength range.

By using components from the LINOS catalog the Q-Sets unify highest precision with a broad area of applications in combination with a flexible handling. The Q-Set components can be used both as kit and individually and allow an easy integration into Microbench setups directly.

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Special Features:

Design facilitates application over a broad wavelength range.

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NEW

Exemplary application for the vertical adjustment set. The base plate is not included in the delivery.

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Q-Set Vertical Adjustment

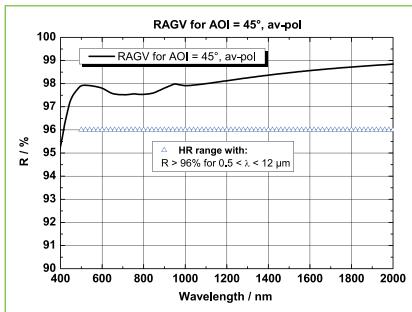
The periscope Q-Set is ideal for applications where the beam height needs to be adapted in a range of 20-180 mm.

Due to the small construction height of the lower beam-deflector cube the system can be easily adapted to the LINOS Microbench level. This enables to integrate a Microbench system in a configuration with a different beam height or vice versa. The maximum beam height is extensible via additional rods.

The mirrors are coated with a high quality silver layer with a reflectivity of more than 96% in the spectral range from 500 nm to 12 μm . Your benefit: They are suitable for applications with broadband light sources, as well as for femtosecond laser pulse applications! The surface-hardened post ensures a very high stability of the structure.

The mirror holder exhibits decoupled adjustment axes allowing a fast and precise alignment process.

The Q-Set is provided as assembly kit:

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Q-Set Variable Attenuator/ Beamsplitter

The Q-Set fulfills two functions, one as variable attenuator where a rotation of the wave plate allows a continuous adjustment of the laser power in a system. Furthermore, it may be used as a variable beam splitter that splits the power between the two orthogonal optical paths.

By its height the Q-Set allows a combination with the LINOS Microbench components. In that case the use of longer rods than the ones included is recommended. In addition, it can be mounted on posts as well as directly onto optical tables.

The zero order wave plates provide an increased wavelength, temperature and incident angle tolerance as wave plates of higher order. For the respective central wavelengths the wave plates are equipped with an anti-reflective coating. The beam splitter cubes are equipped with a broad band anti-reflection coating.

The Q-Set is provided as assembly kit.

Artikel-Nr.	Anz. Qty.	Beschreibung	Description
G065058000	1	Drehfassung	Rotary Support
G061080000	1	Würfel 25	Cube 25
G061207000	4	Stange 20 mm	Rod 20 mm
G061641000	1	Klemmhalter für Prisma 10x10	Clamp holder for prism 10x10
G061253000	1	Basisplatte MB	Mounting Base MB
G335564000 (488 nm) or G335564000 (532 nm) or G335595000 (633 nm) or G335596000 (780 nm)	1	Pol. Strahlteilerwürfel; L=10	Pol. Beamspl. Cube, L=10
G061090000	1	Schraubendreher 1,8 mm	Screw driver 1.8 mm
G061011000	1	Satz Gewindestifte M2,3x3	Set of setscrews M2.3x3 mm
G362701218 (488 nm) or G362701227 (532 nm) or G362701234 (633 nm) or G362701243 (780 nm)	1	Verz.platte D; D=10; L/2; gefasst	Retardation Plate D; D=10; L/2; mounted

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- 10x, NA 0.25, Brennweite 20 mm

Q-Set Spatial Filter

The Q-Set is suited to homogenize the intensity distribution of laser beams. Often dust and dirt particles on optical elements lead to aberrations. Also the generation of the laser light itself may produce an unwanted intensity profile due to cavity properties. These aberrations can be reduced with the aid of the Q-Set Spatial Filter. Here the laser light is focused by an achromatic lens and transmitted through a pinhole which partially removes the inhomogeneities. At the same time the Q-Set can be used to generate a point source.

The pinhole with a diameter of 30 µm is mounted into a x-y-Precision Translation Stage with decoupled adjustment axes. The adjustment screws provide a 0.25 mm spindle pitch and a travel of ±1 mm. The lens can be aligned along the z-axis with an accuracy of 1 µm and almost without any play.

The Q-Set is provided as assembly kit.

XY-translation stage:

- Micrometer screws and fine-adjustment screws with 0.25 mm pitch
- ±1 mm adjustment range
- Decoupled adjustment axis

Z-Fine adjustment:

- Equipped with a micrometer screw with 0.25 mm spindle pitch
- Translation range: 1.5 mm
- Accuracy: 1 µm
- Pinhole: 30 µm diameter

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Q-Set Filter Cascade

The filter cascade provides a stepwise attenuation of light over a huge wavelength and intensity range: the assembly of 5 neutral density filters enables an attenuation over more than 7 orders of magnitude for applications within the Vis and NIR wavelength range. The carriers can be individually flipped in and out of the beam path and therefore allow the desired combination of filters for the respective application. The Q-Set is equipped with neutral density filters with an OD of 0.3; 0.7; 1.3; 2.0 and 3.0. The assembly can be combined with other Microbench setups. Furthermore it allows a mounting directly onto optical tables via the rail which is part of the delivery scope.

The Q-Set is provided as assembly kit.

NEW

Mounting example. (Fibers are not included in the Q-Set.)

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Q-Set Fiber Coupler

The compact design of the Q-Set Fiber Coupler combines a flexible alignment with a high long-term stability. The z-Fine Adjustment unit holding the lens enables a translation of 1 mm with 1 µm resolution and almost without play and tilting. The mounting and alignment of the FSMA fiber coupler is accomplished via a fle.X-plate which is a monolithic element with decoupled adjustment axes in the x and y direction. The respective adjusting range is ±0.25 mm. Adjustment is carried out via tension and pressure screws, which are subsequently secured in position with lock-nuts for excellent long-term stability. The applicability for a broad wavelength range is enabled via the achromatic corrected lens.

The Q-Set is provided as assembly kit.

Fibers are not part of the scope of delivery.

Artikel-Nr. Part No.	Anz. Qty.	Beschreibung	Description
G061209000	4	Stange 75 mm	Rod 75 mm
G061061000	1	Z - Feinverstellung M, Mikrometer	Z-Fine adjustment M micrometer
G038771000	1	Mikroskopobjektiv 10x / 0.25	Microscope objective 10x / 0.25
G067032000	1	fle.X-plate XY Al	fle.X-plate XY Al
G067024000	1	Faseradapter N12,5 FSMA	Fiber optic adapter N12.5 FSMA
G061371000	2	Reiter FLR 40-10	Carrier FLR 40-10
G061352000	1	Flachschiene FLS 40-100	Flat rail FLS 40-100
G061111000	1	Satz Schrauben M2,3x5, 200 Stück	Set of screws M2.3x5, (200pcs.)
G061011000	1	Satz Gewindestifte M2,3x3 (150 Stück)	Set of setscrews M2.3x3 mm (150 pcs.)
G061621000	1	Reduzierring, W0.8" x1/36"	Reducing ring, W0.8" x1/36"
G061090000	1	Schraubendreher 1,8 mm	Screw driver 1.8 mm

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Z-Fine Adjustment:

- No backlash
- High axial precision (almost no tilting)
- Equipped with a micrometer screw with 0.25 mm spindle pitch
- Translation range: 1.5 mm
- Accuracy: 1 µm

Fle.X Plate:

- Monolithic adjustment unit with flexures
- Decoupled adjustment axis
- Position lockable
- Almost no thermal drifting: excellent long-term stability
- The pitch is 0.1 mm per rotation

Q-Set Faserkoppler

Q-Set Fiber Coupler

Produktbezeichnung	Part No.
Item Title	Part No.
Q-Set Fiber Coupler	G060408000

Microbench
nanobench
Tube System C
Positioners
Rail Systems
Mounts and Posts
Mirror Mounts

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Q-Set Spectroscopy Liquids

The Spectroscopy Q-Set for Liquids is designed for use with liquids in standard optic cells with 12.5 mm edge length. Clamping the optic cell to the surface ensures reproducible measurement results.

The modular setup, with components from the Microbench and Flat Rail systems, makes it easy to adapt the structure to any application. See the chapter entitled "Microbench" for compatible components.

The equipment supplied includes a protective cap to prevent light scattering.

Assembled on an FLS Flat Rail 40-100, it can be mounted within a device or on a work surface.

The optic cell and fiber optic cable are not included in the set; please order them separately.

Artikel-Nr. Part No.	Anz. Qty.	Beschreibung	Description
G061081000	1	Würfel 30	Cube 30
G061087000	2	Kollimatorhalter W	Collimator Holder W
G061086000	1	Küvettenhalter	Cuvette Holder
G061078000	2	Verschlusskappe 30	Cap 30
G061352000	1	Schiene FLS 40-100	Rail FLS 40-100
G061374000	1	Reiter FLR 40-40	Carrier FLR 40-40
G120548000	2	SMA Kollimator	SMA Collimator

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The Spectroscopy Q-Set for plates makes it easy to perform transmission and reflection measurements on plates. Mounted with protective plastic screws, it can be used to clamp plates up to 10 mm thick. The collimators can be adjusted to the desired height on the vertical rods, with a Fine Adjustment C30 integrated on one side.

The Q-Set is made up of components of the Microbench and Flat Rail systems and enables individual adaptations and extensions. The foundation of the set is a Flat Rail 40-200, which provides for simple attachment in devices or work surfaces.

The fiber optic cable and glass plate are not included.

Artikel-Nr. Part No.	Anz. Qty.	Beschreibung	Description
G061378000	1	Carrier FLR 40-80	Carrier FLR 40-80
G061374000	2	Carrier FLR 40-40	Carrier FLR 40-40
G061354000	1	Rail FLS 40-200	Rail FLS 40-200
G061085000	3	Mounting Plate FLS	Mounting Plate FLS
G061010000	1	Mounting Plate 25	Mounting Plate 25
G061641000	1	Clamp Holder for Prism 10x10	Clamp Holder for Prism 10x10
G061088000	2	Collimator Holder S	Collimator Holder S
G120548000	2	SMA Collimator	SMA Collimator
G061165000	1	Z-Adjustment M	Z-Adjustment M
G061210000	4	Rail 150 mm	Rod 150 mm
G061207000	4	Rial 20 mm	Rod 20 mm

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Q-Set Retroreflektor

The Q-Set enables the rapid construction of a reflector. The angle between the mirrors can be varied almost arbitrarily and also the distance between input and output beam is variable, whereas even smallest distances are adjustable. An angle fine alignment of the two deflection mirrors is possible via the adjustment screws. The mirror holders are mounted onto a vertically oriented flat rail via carriers. The rail in turn is mounted onto a clamping block attached to a solid steel column and therefore the structure imparts an excellent stability. The distance between the mirrors can be adjusted along the rail to the desired position. The height of the entire system is varied via the clamping block.

In combination with the "Q-Set Prism Stage" the Q-Set Retroreflector is ideal to assemble a prism compressor.

The Q-Set is provided as assembly kit. For assembling three M6x12 screws are required. Mirrors and mounting plate are not included.

Artikel-Nr. Part No.	Anz. Qty.	Beschreibung	Description
G061200000	1	Säule 25 mit Klemmblock	Post 25-200-M with clamping block
G061352000	1	Flachschiene FLS 40-100	Flat Rail FLS 40-100
G061374000	2	Reiter FLR 40-40	Carrier FLR 40-40
G404195000	2	Spiegelhalter MF-25PF	Mirror Mount MF-25PF

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Q-Set Prism Stage

The Q-Set allows a highly accurate and flexible positioning of a prism for different beam heights. The prism table allows an adjustment of the prism in three degrees of freedom about a common pivot point. In addition, a horizontal alignment is possible via the supplied rail (coarse adjustment) and the linear stage (fine adjustment). The mounting column with rod enables adjustment to beam heights 12-21 cm whereas the maximum height can be extended by using longer columns. An ideal area for the Q - set is, for example, together with another identical set in a prism compressor.

The Q-Set is provided as assembly kit. The prism is not included in the delivery. For assembling the following screws are needed : Set screws M6x16 (1x), M2.3x5 (4x); Screw M6x20 (1x), M2.3x3 (4x) M3x6 (4x) M4x8 (4x). These can be found at the end of chapter Microbench and Mounts and Posts.

Artikel-Nr. Part No.	Anz. Qty.	Beschreibung	Description
G061352000	1	Flachschiene FLS 40-100	Flat rail FLS 40-100
G061374000	1	Reiter FLR 40-40	Carrier FLR 40-40
G023041000	1	Säule 14 - 50 - M	Post 14-50-M
G024230000	1	Stift 14x40, 2xM6	Rod 14x40; 2xM6
G061227000	1	Stangenhalter F	Rod Holder F
G061085000	1	Aufnahmeplatte FLS	Mounting Plate FLS
G061207000	2	Stange 20 mm	Rod 20 mm
G085136000	1	Messtisch TB 50-16	Linear stage TB 50-16
G085701000	1	Prismmentisch 40 S	Prism stage 40 S
G061325000	1	Montagewinkel 40	Mounting Bracket

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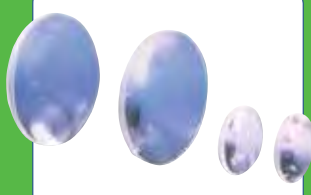
Our quality criteria:

- *Narrowest possible mechanical and optical production tolerances*
- *High-quality coatings*
- *Over 130 years of experience in manufacturing optics*
- *High-quality glass materials from established manufacturers*
- *Available unmounted, or mounted for bench systems*

LINOS Singlets

From convergent lenses and diffuse lenses to best form lenses and aspheres, our extensive selection of simple lenses, or singlets, with various focal lengths and diameters guarantees that you can find the components you need for your application. Qioptiq has over 130 years of experience in the production of precision optical components. Only high-end materials from established manufacturers are used in production, with the narrowest possible production tolerances to ensure the best quality.

Furthermore, the versatile mount system makes LINOS singlets compatible with bench systems both from Qioptiq and from other manufacturers. And, of course, a broad range of economically priced standard anti-reflective coatings is available as well.

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Special Features:

All data (radii, center thicknesses, materials) and designs can also be found in our WinLens optics software packages. Broadband anti-reflective coatings and customer-specific coatings are also available.

Ideal Areas:

Laboratory setups; industrial metrology; laser technology.

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Content

<i>Reflective Curves of Coated Singlets</i>	254
Plano-Convex Lenses	
<i>Plano-Convex Lenses, Unmounted</i>	256
<i>Plano-Convex Lenses, Unmounted (Fused Silica)</i>	258
<i>Plano-Convex Lenses, Mounted</i>	259
<i>Plano-Convex Lenses, Mounted (Fused Silica)</i>	260
Symmetric-Convex Lenses	
<i>Symmetric-Convex Lenses, Unmounted, Dia. 6 - 31.5 mm</i>	262
<i>Symmetric-Convex Lenses, Unmounted (Fused Silica)</i>	263
<i>Symmetric-Convex Lenses, Mounted, Dia. 6 - 22.4 mm</i>	264
<i>Symmetric-Convex Lenses, Mounted, Dia. 25.4 and 31.5 mm</i>	265
<i>Symmetric-Convex Lenses, Mounted (Fused Silica)</i>	266
Plano-Concave Lenses	
<i>Plano-Concave Lenses, Unmounted (N-BK7)</i>	267
<i>Plano-Concave Lenses, Unmounted (Fused Silica)</i>	268
<i>Plano-Concave Lenses, Mounted (N-BK7)</i>	269
<i>Plano-Concave Lenses, Mounted (Fused Silica)</i>	270
Symmetric-Concave Lenses	
<i>Symmetric-Concave Lenses, Unmounted (N-BK7)</i>	271
<i>Symmetric-Concave Lenses, Unmounted (Fused Silica)</i>	272
<i>Symmetric-Concave Lenses, Mounted (N-BK7)</i>	273
<i>Symmetric-Concave Lenses, Mounted (Fused Silica)</i>	274
Aspheric Condenser Lenses	
<i>Aspheric Condenser Lenses, Crown Glass</i>	275
Focus-Variable Lenses from Optotune	
<i>Focus-Variable Lenses from Optotune</i>	276
<i>Electrically Focus Variable Lens EL-10-30</i>	278
<i>Electrically Focus Variable Lens EL-16-40</i>	280
<i>Electrical Lens Driver</i>	281

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Reflective Curves of Coated Singlets

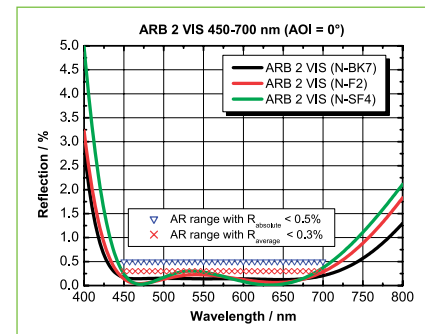
All crown and fused silica singlets of the following pages are also available as coated standard items with rapid delivery.

Broadband Anti-Reflective Coating ARB 2 VIS

- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the visible spectral range*
- *Guaranteed residual reflection $< 0.5\%$ for angles of incidence $0^\circ \leq \text{AOI} \leq 15^\circ$ in the AR*
- *Typical average residual reflection $< 0.3\%$ for angles of incidence $0^\circ \leq \text{AOI} \leq 15^\circ$ in the AR*
- *Damage threshold $H_{\infty} > 10 \text{ J/cm}^2$ at 532 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *AR range typical: 450-700 nm*
- *Design adaptable in the spectral range of $350 \text{ nm} \leq \lambda \leq 950 \text{ nm}$ and for higher angles of incidence*
- *Substrates: N-BK7, fused silica or similar*

Ordering Information:

To order standard lenses with ARB2 VIS on both sides simply change the last three numbers of the order number to 322.

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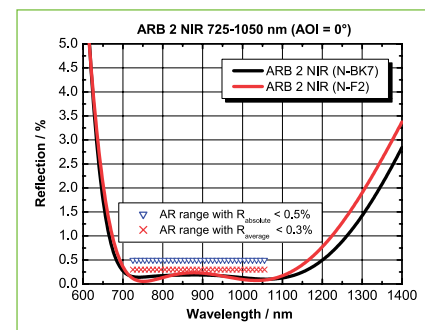
Broadband Anti-Reflective Coating ARB2 NIR

- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the NIR spectral range*
- *Guaranteed residual reflection $< 0.5\%$ for angles of incidence $0^\circ \leq \text{AOI} \leq 15^\circ$ in the AR range*
- *Typical average residual reflection $< 0.3\%$ for angles of incidence $0^\circ \leq \text{AOI} \leq 15^\circ$ in the AR range*
- *Damage threshold: $H_{\infty} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *AR range typical: 725-1050 nm*
- *Design adaptable in the spectral range of $600 \text{ nm} \leq \lambda \leq 1400 \text{ nm}$ and for higher angles of incidence*
- *Substrates: N-BK7, fused silica or similar*

Ordering Information:

To order standard lenses with ARB2 NIR on both sides simply change the last three numbers of the order number to 525.

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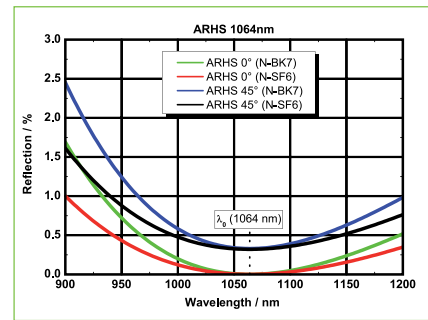
High-power narrowband antireflective coating ARHS

- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for 1064 nm*
- *For Nd:YAG laser applications*

Ordering Information:

To order standard lenses with ARHS 1064 nm on both sides simply change the last three numbers of the order number to 329.

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C*	Optik-Ø (mm)	Sauberkeit ISO 10110	Ø Toleranz (mm)	dm Toleranz (mm)	Zentriergenauigkeit
C*	Ø optics (mm)	Surface Quality ISO 10110	Ø Tolerance (mm)	dm Tolerance (mm)	Centering Accuracy
1	1.5-4	3x0.063	-0.03	±0.1	better than 10'
2	3	3x0.063	-0.09	±0.1	better than 10'
3	5-8	3x0.1	-0.04	±0.1	better than 10'
4	6-8	3x0.063	-0.09	±0.15	better than 4'
5	6-8	3x0.1	-0.09	±0.1	better than 10'
6	10	3x0.1	-0.11	±0.15	better than 4'
7	10	3x0.16	-0.12	±0.15	better than 10'
8	10-18	3x0.1	-0.11	±0.2	better than 4'
9	10-18	3x0.16	-0.12	±0.1	better than 10'
10	12.7	1x0.063	-0.11	±0.2	better than 4'
11	12.7	3x0.16	-0.12	±0.1	better than 10'
12	22.4-31.5	3x0.16	-0.13	±0.2	better than 4'
13	22.4-31.5	3x0.25	-0.15	±0.2	better than 10'
14	25.4-31.5	3x0.25	-0.07	±0.2	better than 3'
15	40-50	5x0.25	-0.19	±0.2	better than 10'
16	50	3x0.25	-0.16	±0.2	better than 4'

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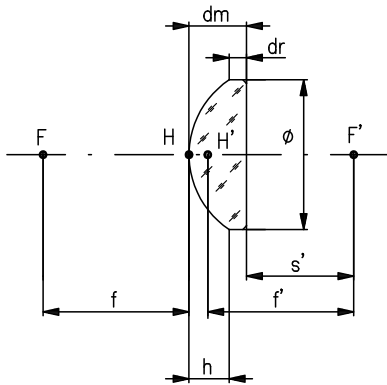
Our lenses are also available with the following standard, economically priced anti-reflective coatings:

- ARB 2 VIS (450-700 nm), order code -322
- ARB 2 NIR (725-1050 nm), order code -525
- ARHS 1064 (Nd:YAG center wavelength at 1064 nm), order code -329

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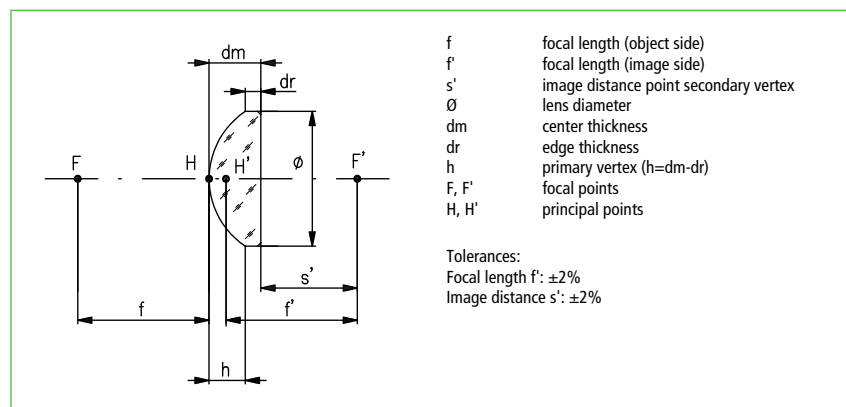
Plano-Convex Lenses, Unmounted (Fused Silica)

- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$

Plankonvexlinsen, ungefasst, aus Quarzglas Plano-convex lenses, unmounted, fabricated from fused silica

Brennweite (mm)	Ø (mm)	d_m (mm)	d_r (mm)	$s'_{248\text{ nm}}$ (mm)	$f'_{248\text{ nm}}$ (mm)	C^*	Part No.
Focal length (mm)	Ø (mm)	d_m (mm)	d_r (mm)	$s'_{248\text{ nm}}$ (mm)	$f'_{248\text{ nm}}$ (mm)	C^*	Part No.
12.5	8	2.4	1	10.91	12.5	4	G312255000
15	10	2.9	1	12.83	14.75	8	G312256000
20	12.5	3.5	1.3	17.35	19.67	8	G312257000
25	12.5	3	1.3	22.78	24.77	8	G312258000
30	18	4.4	1.5	27.59	30.51	8	G312259000
40	22.4	5	1.5	35.38	38.69	12	G312286000
50	22.4	4.1	1.5	46.7	49.42	12	G312300000
60	22.4	3.7	1.6	57.13	59.58	12	G312301000
80	22.4	3.1	1.5	77.4	79.45	12	G312302000
100	22.4	2.8	1.5	97.45	99.31	12	G312303000
200	22.4	2.1	1.5	195.33	196.73	12	G312304000
40	31.5	9.5	2	33.65	39.95	12	G312411000
50	31.5	7.5	2.1	45.14	50.11	12	G312412000
60	31.5	6.4	2	55.75	60	12	G312413000
80	31.5	5.2	2	76.55	80	12	G312414000
100	31.5	4.5	2	96.99	99.97	12	G312415000
140	31.5	3.3	1.5	136.09	138.21	12	G312305000
200	31.5	3.3	2.1	198.03	200.22	12	G312416000
300	31.5	3	2.2	298.66	300.65	12	G312417000
400	31.5	2.1	1.5	390.96	392.35	12	G312306000
100	50	8.5	1.8	93.67	99.31	16	G312242000
150	50	6.1	1.8	144.54	148.59	16	G312243000
200	50	5.1	1.9	193.34	196.73	16	G312244000
250	50	4.4	1.9	244.74	247.66	16	G312245000

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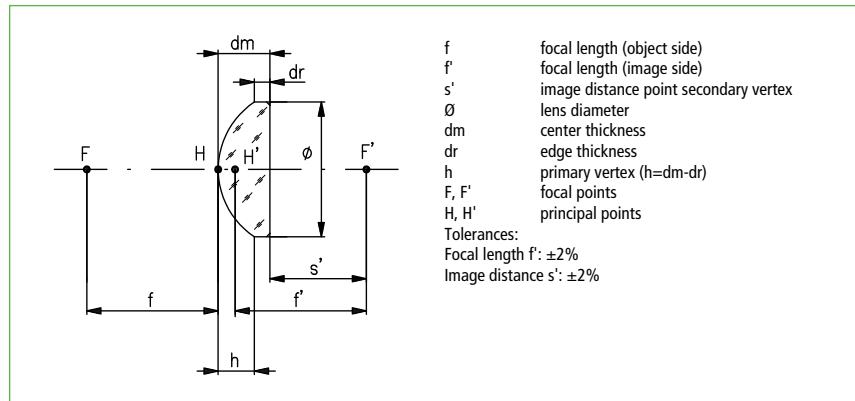
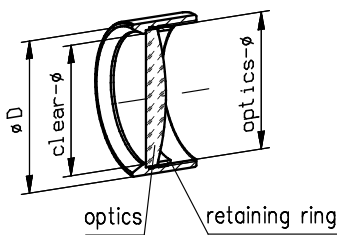
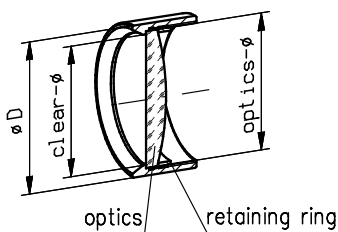
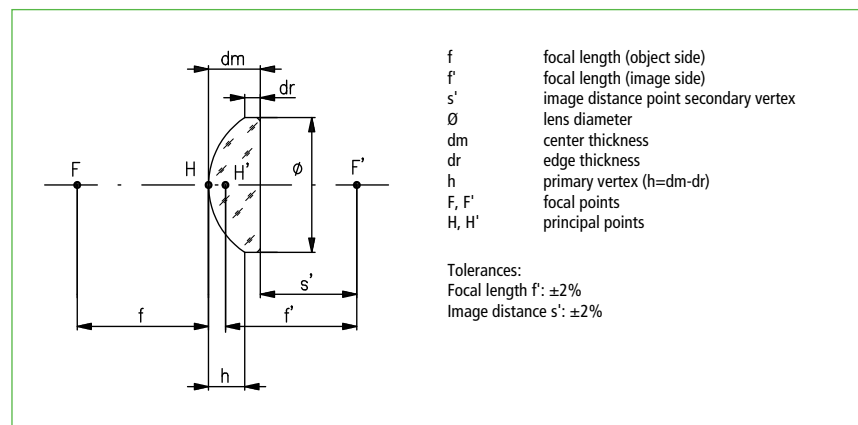
All relevant specifications of the lenses, such as materials, radii and center thicknesses, are included in the database of the our WinLens optical design software. A freeware version of this program is available for download at winlens.de.

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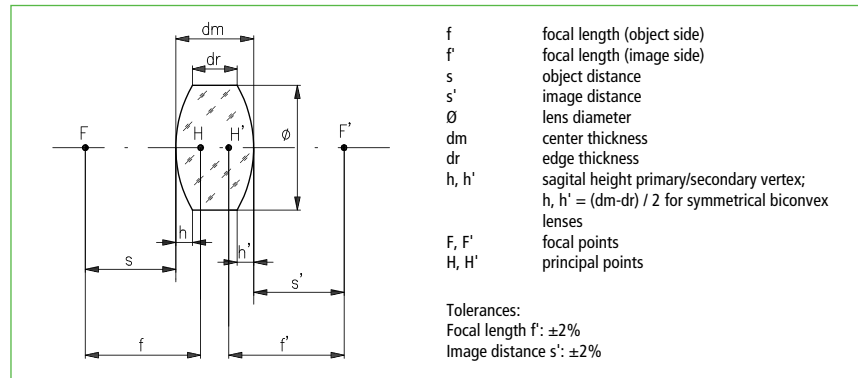
Symmetric-Convex Lenses, Unmounted, Dia. 6 - 31.5 mm

- Precision designed for 1:1 imaging
- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$
- Two equal radii of curvature

Bikonvexlinsen, ungefasst, Durchmesser 6 - 31.5 mm Symmetric-convex lenses, unmounted, dia. 6 - 31.5 mm


Brennweite (mm)	\emptyset (mm)	Material	dm (mm)	dr (mm)	$s, s'_{588\text{ nm}}$ (mm)	$f'_{546\text{ nm}}$ (mm)	C^*	Part No.
Focal length (mm)	\emptyset (mm)	Material	dm (mm)	dr (mm)	$s, s'_{588\text{ nm}}$ (mm)	$f'_{546\text{ nm}}$ (mm)	C^*	Part No.
5	6	N-BaK4	2.9	0.9	3.95	4.97	5	G311328000
7	8	N-BaK4	3.4	1.1	6	7.14	5	G311329000
10	10	N-BaK4	3.5	1	8.78	9.93	9	G311300000
12.5	12.5	N-BK7	4.5	1	10.79	12.33	9	G311377000
15	12.7	N-BK7	4	1.1	13.58	14.91	9	G311381000
20	12.7	N-BK7	3	0.9	18.94	19.88	9	G311382000
30	12.7	N-BK7	3	1.6	28.8	29.7	9	G311383000
50	12.7	N-BK7	3	2.2	48.92	49.74	9	G311384000
150	12.7	N-BK7	3	2.7	148.5	148.9	9	G311387000
16	18	N-BK7	7	1.1	13.43	15.88	9	G311378000
20	22.4	N-BK7	8.8	1.6	17.05	20.13	13	G311338000
25	22.4	N-BK7	7	1.6	22.4	24.74	13	G311308000
30	22.4	N-BK7	6.3	2	28.01	30.06	13	G311310000
40	22.4	N-BK7	5	2	39.66	41.2	13	G311314000
50	22.4	N-BK7	4.5	2.1	49.88	51.19	13	G311315000
60	22.4	N-BK7	4	2	59.94	61.05	13	G311316000
80	22.4	N-BK7	3.5	2	80.24	81.1	13	G311317000
100	22.4	N-BK7	3.5	2.3	100.44	101.22	13	G311318000
150	22.4	N-BK7	3	2.2	150.65	151.08	13	G311319000
200	22.4	N-BK7	3	2.4	202.51	202.75	13	G311320000
30	25.4	N-BK7	7	1.4	27.89	30.18	14	G311321000
50	25.4	N-BK7	7	3.8	48.24	50.42	14	G311322000
80	25.4	N-BK7	7	5	77.36	79.41	14	G311323000
100	25.4	N-BK7	7	5.4	98.41	100.4	14	G311324000
150	25.4	N-BK7	7	6	147.8	149.6	14	G311325000
30	31.5	N-SK2	9.5	2	27.03	30.02	14	G311331000
40	31.5	N-BK7	8.3	2.1	39.09	41.77	14	G311339000
50	31.5	N-BK7	7	2.1	49.45	51.62	14	G311351000
70	31.5	N-BK7	5.5	2.1	69.57	71.15	14	G311357000
100	31.5	N-BK7	4.8	2.4	98.41	100	14	G311333000
130	31.5	N-BK7	4	2.2	129.3	130.14	14	G311335000

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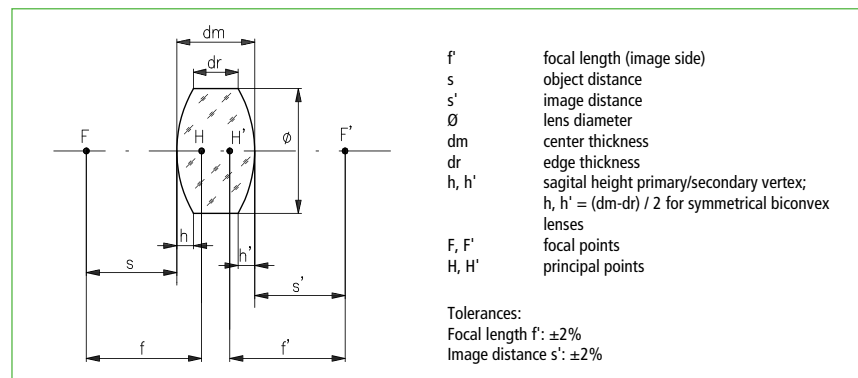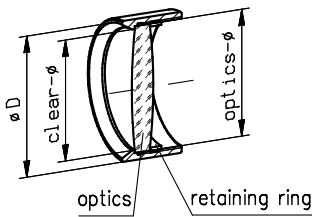
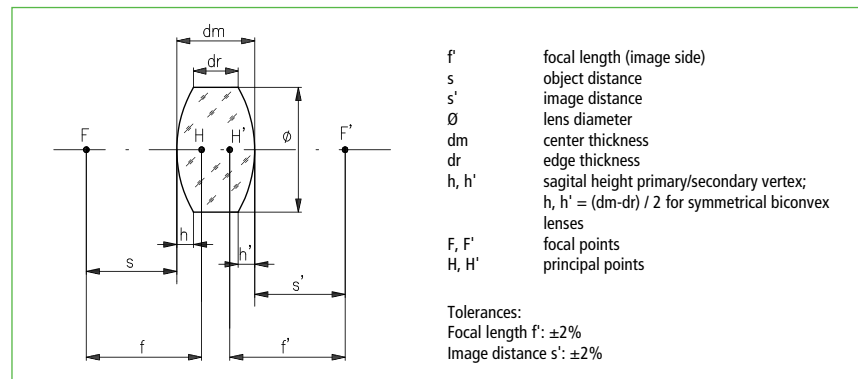
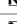
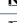
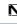
Symmetric-Convex Lenses, Mounted, Dia. 6 - 22.4 mm


- Lens mounts fit to Microbench  or Nanobench 
- Precision designed for 1:1 imaging
- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$
- Two equal radii of curvature

Bikonvexlinsen, gefasst, Durchmesser 6 - 22.4 mm Symmetric-convex lenses, mounted, dia. 6 - 22.4 mm

Fassung	Brennweite (mm)	Optik- ϕ (mm)	Material	Außen- ϕ D (mm)	Freier- ϕ (mm)	dm (mm)	dr (mm)	$s, s'_{588\text{ nm}}$ (mm)	$f'_{546\text{ nm}}$ (mm)	ϕ^*	Part No.
Mount	Focal length (mm)	ϕ optics (mm)	Material	Outer- ϕ D (mm)	Clear ϕ (mm)	dm (mm)	dr (mm)	$s, s'_{588\text{ nm}}$ (mm)	$f'_{546\text{ nm}}$ (mm)	ϕ^*	Part No.
	5	6	N-BaK4	16	5	2.9	0.9	3.95	4.97	5	G052201000
	7	8	N-BaK4	16	7	3.4	1.1	6	7.14	5	G052202000
	10	10	N-BaK4	16	9	3.5	1	8.78	9.93	9	G052203000
	12.5	12.5	N-BK7	16	11.5	4.5	1	10.79	12.33	9	G052204000
	15	12.7	N-BK7	16	11.5	4	1.1	13.58	14.91	9	G052205000
	20	12.7	N-BK7	16	11.5	3	0.9	18.94	19.88	9	G052206000
	30	12.7	N-BK7	16	11.5	3	1.6	28.8	29.7	9	G052207000
	50	12.7	N-BK7	16	11.5	3	2.2	48.92	49.74	9	G052208000
	100	12.7	N-BK7	16	11.5	3	2.6	99.08	99.7	9	G052210000
	150	12.7	N-BK7	16	11.5	3	2.7	148.5	148.9	9	G052211000
	16	18	N-BK7	25	17	7	1.1	13.43	15.88	9	G063033000
	20	22.4	N-BK7	25	21.4	8.8	1.6	17.05	20.13	13	G063020000
	25	22.4	N-BK7	25	21.4	7	1.6	22.4	24.74	13	G063021000
	30	22.4	N-BK7	25	21.4	6.3	2	28.01	30.06	13	G063022000
	40	22.4	N-BK7	25	21.4	5	2	39.66	41.2	13	G063023000
	50	22.4	N-BK7	25	21.4	4.5	2.1	49.88	51.19	13	G063024000
	60	22.4	N-BK7	25	21.4	4	2	59.94	61.05	13	G063025000
	80	22.4	N-BK7	25	21.4	3.5	2	80.24	81.1	13	G063026000
	100	22.4	N-BK7	25	21.4	3.5	2.3	100.44	101.22	13	G063027000
	150	22.4	N-BK7	25	21.4	3	2.2	150.65	151.08	13	G063028000
	200	22.4	N-BK7	25	21.4	3	2.4	202.51	202.75	13	G063029000

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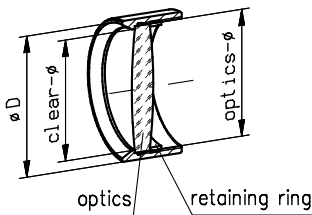
Symmetric-Convex Lenses, Mounted, Dia. 25.4 and 31.5 mm

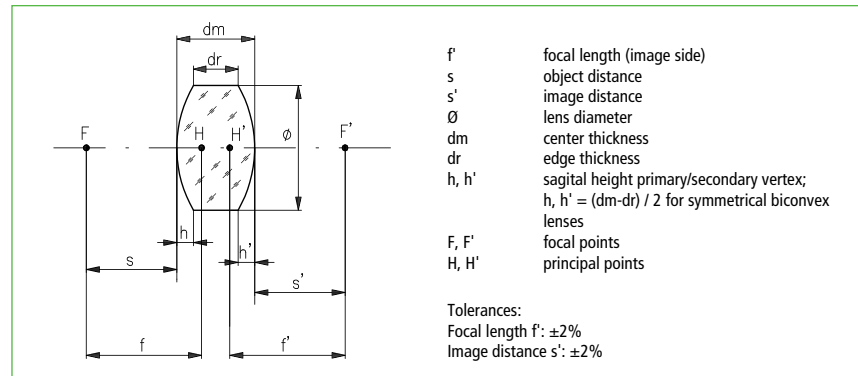- Lens mounts fit to Microbench ☒
- Precision designed for 1:1 imaging
- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$
- Two equal radii of curvature

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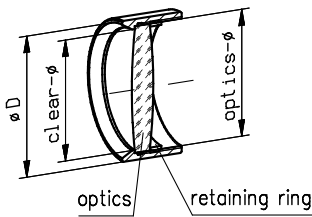
For specifications of the mounts, such as dimensions, tolerances, etc., please look at chapter "Mounts and Posts", section "Mounts for Microbench and Nanobench".

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
Symmetric-Convex Lenses, Mounted (Fused Silica)

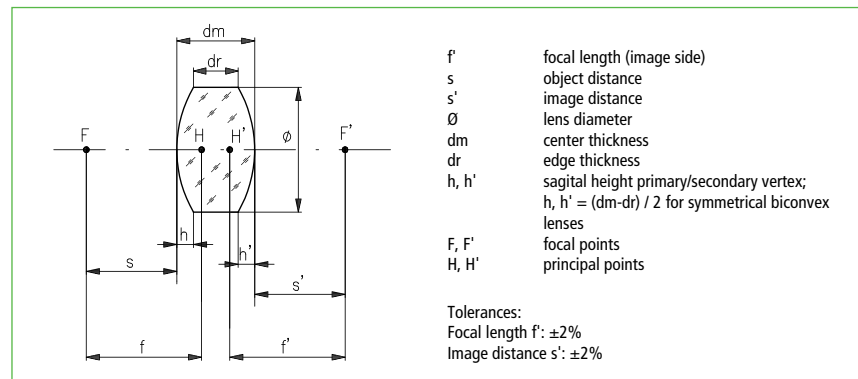- Lenses mounts fit to Microbench or Nanobench
- Precision designed for 1:1 imaging
- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$
- Two equal radii of curvature

Bikonvexlinsen, gefasst, aus Quarzglas Symmetric-convex lenses, mounted, fabricated from fused silica

Fassung	Brennweite (mm)	Optik- \emptyset (mm)	Außen- \emptyset D (mm)	Freier- \emptyset (mm)	dm (mm)	dr (mm)	$S, S'_{248\text{ nm}}$ (mm)	$f'_{193\text{ nm}}$ (mm)	$f'_{248\text{ nm}}$ (mm)	C^*	Part No.
Mount	Focal length (mm)	\emptyset optics (mm)	Outer- \emptyset D (mm)	Clear \emptyset (mm)	dm (mm)	dr (mm)	$S, S'_{248\text{ nm}}$ (mm)	$f'_{193\text{ nm}}$ (mm)	$f'_{248\text{ nm}}$ (mm)	C^*	Part No.
	5	6	16	5	3	1	4.38	5.01	5.48	4	G052212000
	10	10	16	9	4	1.2	8.71	9.24	10.13	8	G052213000
	12.5	12.5	16	11.5	4.7	1	10.52	11.11	12.2	8	G052214000
	16	12.5	16	11.5	3.8	1.2	14.82	14.67	16.13	8	G052215000
	20	18	25	17	5.8	1.4	18.33	18.52	20.36	8	G063316000
	25	22.4	25	21.4	7	1.6	22.79	22.95	25.23	12	G063030000
	30	22.4	25	21.4	6	1.5	27.5	26.87	29.56	12	G063317000
	40	22.4	25	21.4	4.9	1.7	38.05	36.05	39.71	12	G063318000
	50	22.4	25	21.4	4.2	1.6	47.89	44.74	49.3	12	G063319000
	60	22.4	25	21.4	3.7	1.6	58.82	54.49	60.06	12	G063320000
	80	22.4	25	21.4	3.3	1.7	76.47	70.36	77.57	12	G063031000
	100	22.4	25	21.4	2.9	1.6	97.88	89.65	98.85	12	G063321000
	150	22.4	25	21.4	2.4	1.6	146.77	133.82	147.57	12	G063322000
	200	22.4	25	21.4	2.2	1.6	195.9	178.3	196.63	12	G063323000
	30	31.5	35	30	11.5	2	25.97	27.39	30.06	12	G063161000
	40	31.5	35	30	8.6	2	37.1	36.43	40.06	12	G063162000
	50	31.5	35	30	7.3	2.2	47.52	45.42	50	12	G063163000
	60	31.5	35	30	6.2	2	57.95	54.51	60.04	12	G063164000
	80	31.5	35	30	5.1	2	78.37	72.68	80.08	12	G063165000
	100	31.5	35	30	4.5	2	98.64	90.86	100.14	12	G063166000
	120	31.5	35	30	4.1	2.1	118.72	108.95	120.09	12	G063168000

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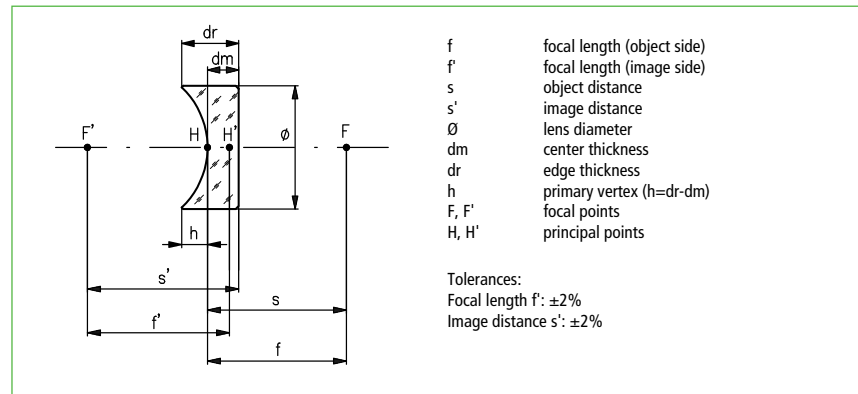
Plano-Concave Lenses, Unmounted (N-BK7)

- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$

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All relevant specifications of the lenses, such as materials, radii and center thicknesses, are included in the database of our LINOS WinLens optical design software. A freeware version of this program is available for download at winlens.de.

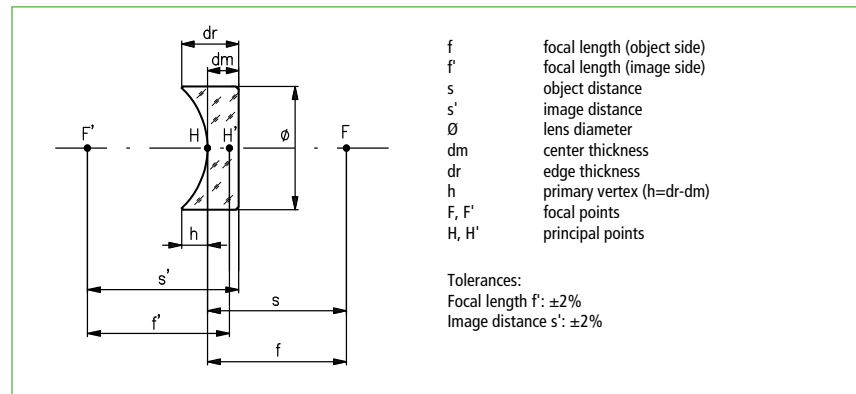
- Informationen zu Laserspiegelschichten finden Sie im Kapitel "Dünne Schichten"
- For information on laser mirror coatings, please refer to chapter "Thin Film Coatings"

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Plano-Concave Lenses, Unmounted (Fused Silica)

- Substrate for dielectric or metallic coatings
- Suitable for DUV applications
- Uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$
- Extreme low scattering losses
- High surface quality
- Surface quality: $\lambda/10$ for concave surfaces, $\lambda/4$ for flat surfaces (P-V).

Plankonkavlinen, ungefasst (aus Quarzglas) Plano-concave lenses, unmounted (fabricated from fused silica)

Brennweite (mm)	\varnothing (mm)	dm (mm)	dr (mm)	$s'_{248\text{ nm}}$ (mm)	$s_{248\text{ nm}}$ (mm)	$f'_{193\text{ nm}}$ (mm)	$f'_{248\text{ nm}}$ (mm)	C^*	Part No.
Focal length (mm)	\varnothing (mm)	dm (mm)	dr (mm)	$s'_{248\text{ nm}}$ (mm)	$s_{248\text{ nm}}$ (mm)	$f'_{193\text{ nm}}$ (mm)	$f'_{248\text{ nm}}$ (mm)	C^*	Part No.
-10	6	1	2	-10.52	9.36	-8.94	-9.85	4	G314419000
-16	10	1	2.7	-16.74	15.28	-14.58	-16.08	6	G314420000
-20	12.7	4	6	-22.32	19.66	-17.83	-19.66	10	G340160000
-40	22.4	1.5	4.9	-40.51	37.55	-35.84	-39.52	12	G314421000
-50	22.4	1.5	4.1	-51.1	47.62	-45.44	-50.11	12	G314422000
-100	22.4	1.5	2.8	-100.26	94.33	-90.02	-99.27	12	G314423000
-150	22.4	1.5	2.3	-149.51	141.14	-134.69	-148.52	12	G314424000

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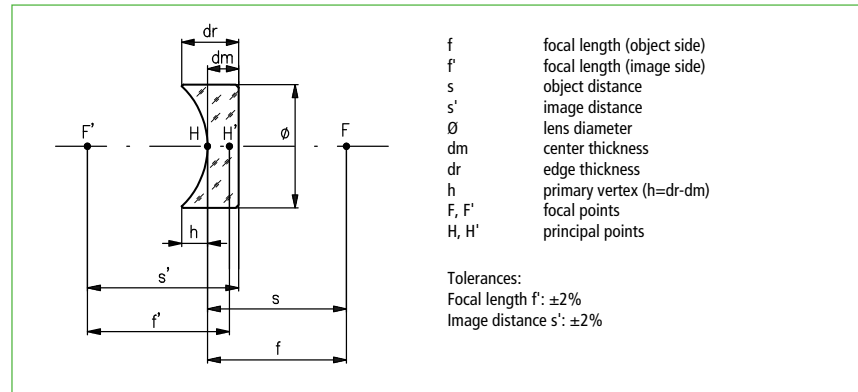
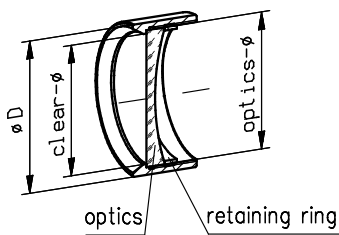
Plano-Concave Lenses, Mounted (N-BK7)

- Mounted lenses fit to Microbench ☒ or Nanobench ☒
- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$

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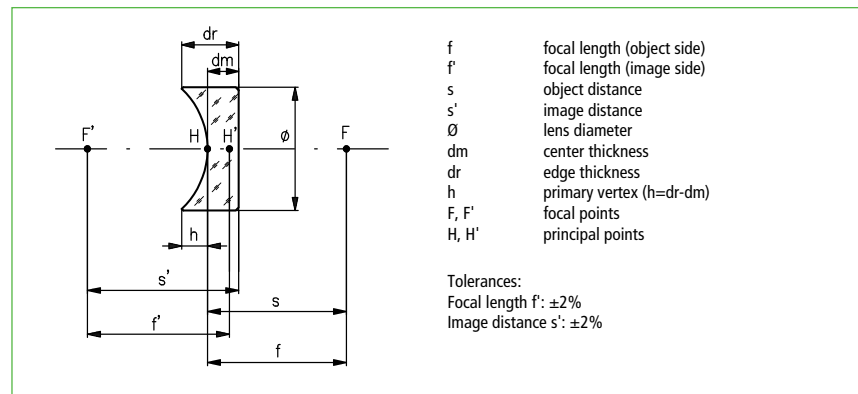
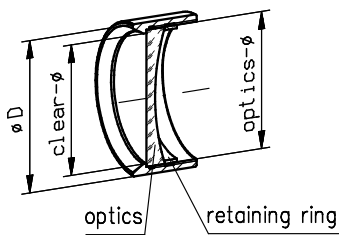
Plano-Concave Lenses, Mounted (Fused Silica)


- Mounted lenses fit to Microbench ☒ or Nanobench ☒
- Suitable for DUV applications
- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$

Plankonkavlinen, gefasst aus Quarzglas Plano-concave lenses, mounted (fabricated from fused silica)

Fassung	Brennweite (mm)	Optik- ϕ (mm)	Außen- ϕD (mm)	Freier- ϕ (mm)	dm (mm)	dr (mm)	$s'_{248\text{ nm}}$ (mm)	$f'_{248\text{ nm}}$ (mm)	C^*	Part No.
Mount	Focal length (mm)	ϕ optics (mm)	Outer- ϕD (mm)	Clear ϕ (mm)	dm (mm)	dr (mm)	$s'_{248\text{ nm}}$ (mm)	$f'_{248\text{ nm}}$ (mm)	C^*	Part No.
☒	-10	6	16	5	1	2	-10.52	-9.85	4	G052313000
☒	-16	10	16	9	1	2.7	-16.74	-16.08	8	G052314000
☒	-20	12.7	16	11.5	4	6	-22.32	-19.66	8	G052315000
☒	-200	12.7	16	11.5	5.8	6	-200.48	-196.64	8	G052318000
☒	-40	22.4	25	21.4	1.5	4.9	-40.51	-39.52	12	G063352000
☒	-50	22.4	25	21.4	1.5	4.1	-51.1	-50.11	12	G063353000
☒	-100	22.4	25	21.4	1.5	2.8	-100.26	-99.27	12	G063354000
☒	-150	22.4	25	21.4	1.5	2.3	-149.51	-148.52	12	G063355000

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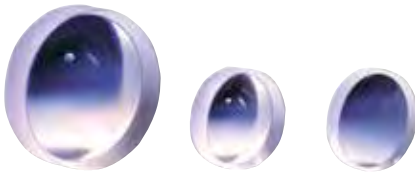
Symmetric-Concave Lenses, Unmounted (N-BK7)

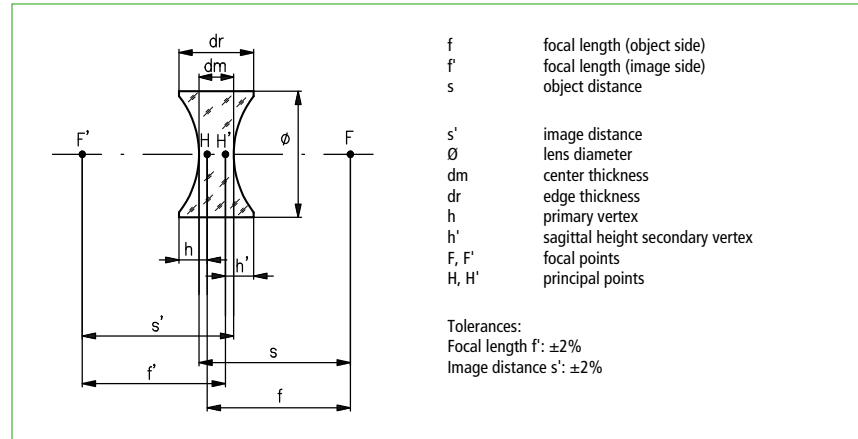- Two equal radii of curvature
- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$

Bikonkavlinen, ungefasst (N-BK7), symmetrisch Symmetric-concave lenses, unmounted (N-BK7)


Brennweite (mm)	\varnothing (mm)	dm (mm)	dr (mm)	h, h' (mm)	$s, s'_{588\text{ nm}}$ (mm)	$f'_{546\text{ nm}}$ (mm)	C^*	Part No.
Focal length (mm)	\varnothing (mm)	dm (mm)	dr (mm)	h, h' (mm)	$s, s'_{588\text{ nm}}$ (mm)	$f'_{546\text{ nm}}$ (mm)	C^*	Part No.
-10	10	1	3.5	1.24	-10.48	-10.12	7	G313210000
-20	12.7	1.5	3.5	0.98	-20.55	-19.98	11	G313373000
-50	12.7	1.5	2.3	0.39	-50.88	-50.2	11	G313374000
-100	12.7	1.5	1.9	0.19	-101.3	-100.4	11	G313375000
-20	22.4	1.5	7.5	3.27	-20.4	-19.84	13	G313205000
-30	22.4	1.5	5.6	2.07	-30.62	-30.02	13	G313206000
-40	22.4	1.5	4.5	1.53	-40.75	-40.11	13	G313214000
-50	22.4	1.5	3.9	1.21	-50.87	-50.2	13	G313215000
-100	22.4	1.5	2.7	0.6	-101.26	-100.4	13	G313216000
-150	22.4	1.5	2.3	0.4	-151.39	-150.34	13	G313217000

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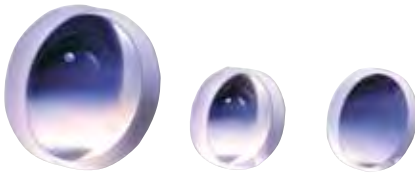
Symmetric-Concave Lenses, Unmounted (Fused Silica)

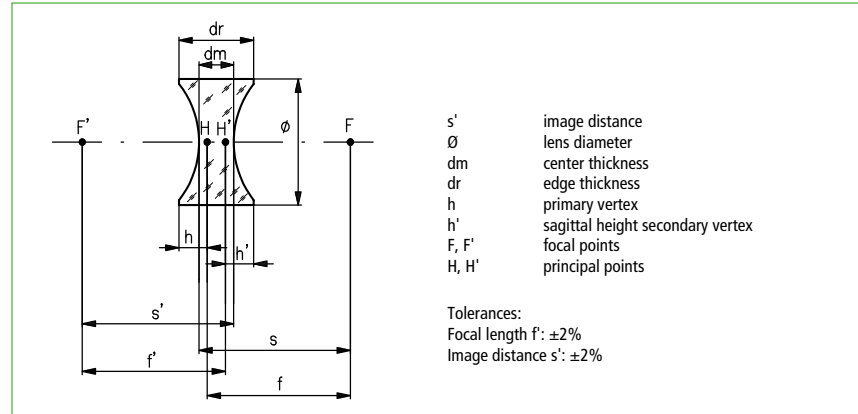- Suitable for DUV-applications
- Two equal radii of curvature
- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$

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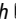
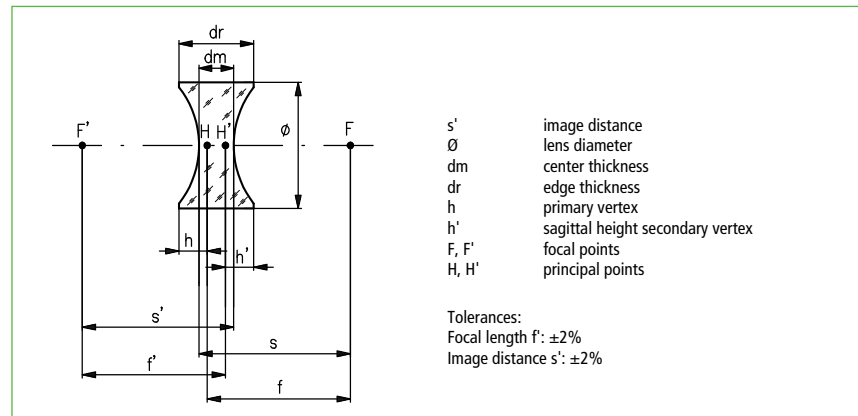
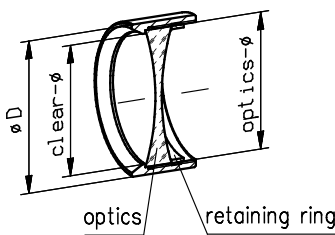
Symmetric-Concave Lenses, Mounted (N-BK7)

- Two equal radii of curvature
- Mounted lenses fit to Microbench  or Nanobench 
- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$

Bikonkavlinen, gefasst (N-BK7), symmetrisch Symmetric-concave lenses, mounted (N-BK7)

Fassung	Brennweite (mm)	Optik- \varnothing (mm)	Außen- $\varnothing D$ (mm)	Freier- \varnothing (mm)	dm (mm)	dr (mm)	h, h' (mm)	$s, s'_{588\text{ nm}}$ (mm)	$f'_{546\text{ nm}}$ (mm)	C^*	Part No.
Mount	Focal length (mm)	\varnothing optics (mm)	Outer- $\varnothing D$ (mm)	Clear \varnothing (mm)	dm (mm)	dr (mm)	h, h' (mm)	$s, s'_{588\text{ nm}}$ (mm)	$f'_{546\text{ nm}}$ (mm)	C^*	Part No.
	-10	10	16	9	1	3.5	1.24	-10.48	-10.12	7	G052401000
	-20	12.7	16	11.5	1.5	3.5	0.98	-20.55	-19.98	11	G052402000
	-50	12.7	16	11.5	1.5	2.3	0.39	-50.88	-50.2	11	G052403000
	-100	12.7	16	11.5	1.5	1.9	0.19	-101.3	-100.4	11	G052404000
	-20	22.4	25	21.4	1.5	7.2	3.27	-20.4	-19.84	13	G063058000
	-30	22.4	25	21.4	1.5	5.6	2.07	-30.62	-30.02	13	G063059000
	-40	22.4	25	21.4	1.5	4.5	1.53	-40.75	-40.11	13	G063060000
	-50	22.4	25	21.4	1.5	3.9	1.21	-50.87	-50.2	13	G063061000
	-100	22.4	25	21.4	1.5	2.7	0.6	-101.26	-100.4	13	G063062000
	-150	22.4	25	21.4	1.5	2.3	0.4	-151.39	-150.34	13	G063063000

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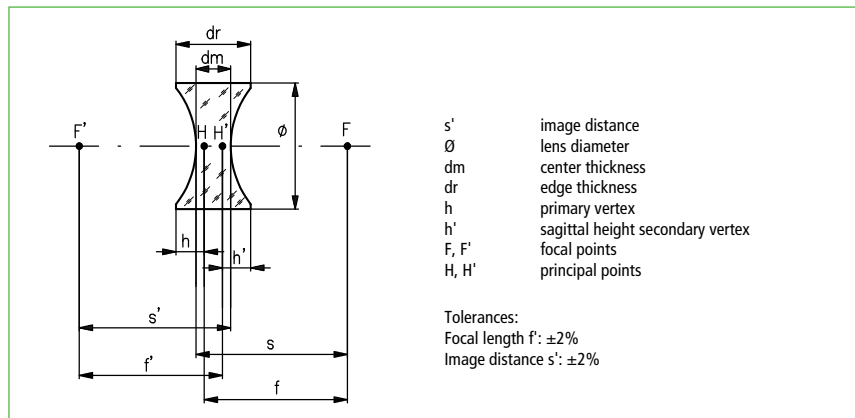
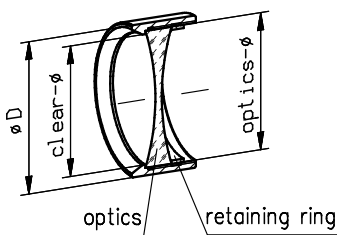
Symmetric-Concave Lenses, Mounted (Fused Silica)

- Suitable for DUV applications
- Two equal radii of curvature
- Mounted lenses fit to Microbench
- Surfaces uncoated
- Tolerances: Focal length f' : $\pm 2\%$, Image distance s' : $\pm 2\%$

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
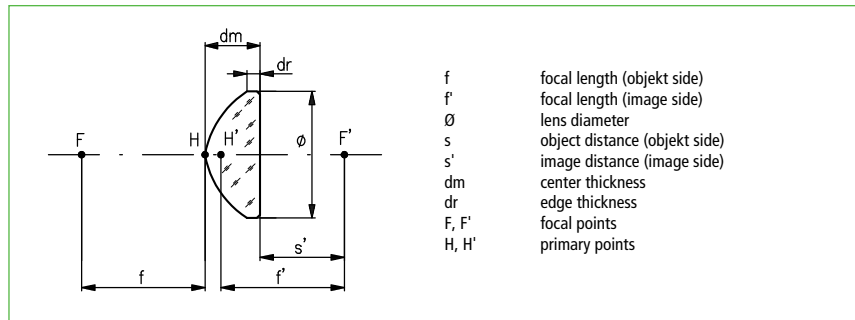
Aspheric Condenser Lenses, Crown glass

- *Plano-convex lenses with minimal spherical aberration*
- *Very high aperture ratios*
- *Primarily for use in illumination systems*
- *Mounts fit to Microbench ☒*
- *Fabricated from crown glass*
- *Surfaces uncoated*

Asphärische Kondensorenlinen, ungefasst (Kronglas) Aspheric condenser lenses, unmounted (crown glass)

Brennweite (mm)	Ø (mm)	dm (mm)	dr (mm)	Ø Toleranz (mm)	$s'_{588\text{ nm}}$ (mm)	$f'_{546\text{ nm}}$ (mm)	$f'_{588\text{ nm}}$ (mm)	Part No.
15	18	7.4	2	-0.13	10.7	15.7	15.8	G317703000
18	22.4	10.2	2.5	-0.13	11.8	18.7	18.8	G317704000
27	31.5	12.3	1.8	-0.16	19.5	27.8	27.9	G317707000
40	50	19.4	2.3	-0.3	28.6	41.8	42	G317708000

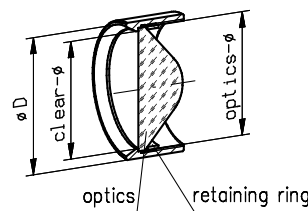Asphärische Kondensorenlinen, gefasst (Kronglas) Aspheric condenser lenses, mounted (crown glass)

Fassung	Brennweite (mm)	Ø (mm)	Fassung, Außen-Ø D (mm)	dm (mm)	dr (mm)	Ø Toleranz (mm)	$s'_{588\text{ nm}}$ (mm)	$f'_{546\text{ nm}}$ (mm)	$f'_{588\text{ nm}}$ (mm)	Part No.
☒	15	18	25 mm, clear Ø 17 mm	7.4	2	-0.13	10.7	15.7	15.8	G063095000
☒	18	22.4	25 mm, clear Ø 21.4 mm	10.2	2.5	-0.13	11.8	18.7	18.8	G063097000
☒	27	31.5	35 mm, clear Ø 30 mm	12.3	1.8	-0.16	19.5	27.8	27.9	G063098000

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Focus-Variable Lenses from Optotune

Functionality and characteristics

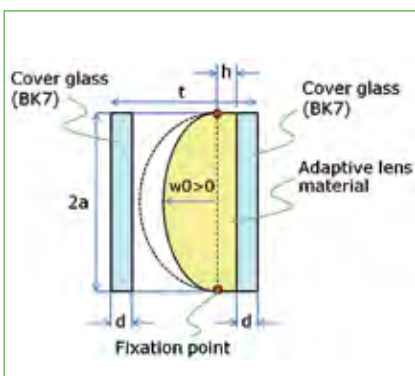
The focus-variable lenses from Optotune are manufactured from elastic polymers and have unique properties: the lens shape can be changed from convex to concave mechanically or electronically. A clear aperture from 1 to 55 mm is possible.

Both lens types work with the same principle: two chambers are separated through a thin membrane, one chamber is filled with a liquid and the other with air. The pressure difference between the two chambers defines the deflection and with that the radius of the lens. The pressure difference is being controlled by using a thread ring to push a ring-shaped actuator down onto the chamber. The changes are reversible; there is no hysteresis.

Advantages

The variable focal length of these innovative spherical lenses enables the design of very compact optical systems. A change in the radius of curvature of just a few micrometers can correspond to an axial shift in the centimeter range. The typical switching frequency of the electronically adjustable lens is of the order of about 100 Hz.

Fewer lenses will be needed for an optical design; an axial movement of the components is less complex or even unnecessary. No need for complex mechanics or motors anymore. This means: fewer moving components and thus a lighter, more rugged, fully closed and dust-free design.

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- a: lens radius
- w0: central bending (depends on current)
- h: constant zone of the lense material = 2.45 ± 0.3 mm (EL-10-30)
- d: thickness of the cover glass (BK7) = 0.5 mm
- t: overall thickness of the lense = 8.5 mm ± 0.3 mm (EL-10-30)
- red: fixing of the membrane
- yellow: adaptive lens material (Optotune)

Optical design of the electrically focus tunable lense EL-10-30

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Field of Application

These lenses are ideally suited to be used both in compact imaging systems, laser material processing and in illumination systems, such as LED-based systems.

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Compact Zoom Optic

A change of the radius of curvature of a lens has a greater visual effect than is common practice with adjusting lenses. This enables the development of more compact optical systems. This way it is possible to save up to 50% of the overall length of e.g. an optical 3-times zoom lens.

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Fast Autofocus

Applications in the Machine Vision field often demand very fast focusing onto different working distances. This most often requires that the radius of one single lens is being adjusted – which is the ideal application area of the focus-variable lens.

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About the Manufacturer

The Optotune AG is a spin-off of the ETH Zurich. It is active in research, development and production of adaptive optical systems. The core competence is processing the used materials, especially polymers, and to develop more compact optical systems.

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Advantages of focus-variable lenses from Optotune:

- *Compact design*
- *Fast response time*
- *Robustness (no complex mechanics)*
- *High switching frequency (frictionless)*

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Electrically Focus Variable Lens EL-10-30

Compact optics and short setting time: EL-10-30 allows for that due to its electrically tunable focal range. Within 10 ms, the radius of this plano-convex lens can be changed by applying a voltage between 0 and 5 V. The lens can be operated with commercial power supplies allowing for current control. Power consumption is less than 2 W. For each lens a test chart displaying the dependency between focal length and current can be provided.



The EL-10-30 comes in two different housings: Compact and C-mount. While the compact housing is preferably mounted in a 30 mm ring, the C-mount housing has threads on either side. Furthermore, the C-mount version accommodates space to insert an offset lens of 12 or 12.7mm diameter. The working principle of both versions is identical. C-mount versions marked with a "C" in the name.

In the download area, a DLL is provided for embedding the lens in ZEMAX.

- Aperture 10 mm
- Lifecycles (10-90% sinusoidal): >1.000.000.000
- Response time (10%-90% step): < 2.5 ms
- Settling time < 15 ms
- Polarization preserving
- Transmission > 90 % @ 633 nm
- AR coated for 400-700 nm or 700-1100 nm
- Abbe number: 31 (HR) and 100 (LD)
- HR:
 - n_F (486.1 nm) = 1.572,
 - n_D (589.3 nm) = 1.559,
 - n_C (656.3 nm) = 1.554
- LD:
 - n_F (486.1 nm) = 1.302,
 - n_D (589.3 nm) = 1.300,
 - n_C (656.3 nm) = 1.299
- Damage threshold 10 kW/cm² (@1064 nm)
- Fits to Microbench with holder G06 1042 000 or G06 1225 000

Elektrisch Fokus-Variable Linse EL-10-30 Electrically Focus Variable Lens EL-10-30


Produktbezeichnung	Brennweitenbereich	Part No.
Item Title	Focal Range	Part No.
EL-10-30-TC-VIS-12D	+50 to +120	G399553100
EL-10-30-TC-NIR-12D	+50 to +120	G399554100
EL-10-30-Ci-VIS-LD	+80 to +200	G399569000
EL-10-30-Ci-NIR-LD	+80 to +200	G399571000
EL-10-30-Ci-VIS-LD-MV	-600 to +170	G399572000

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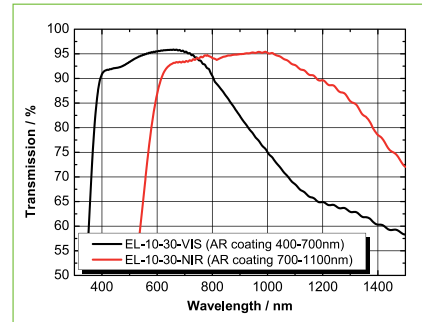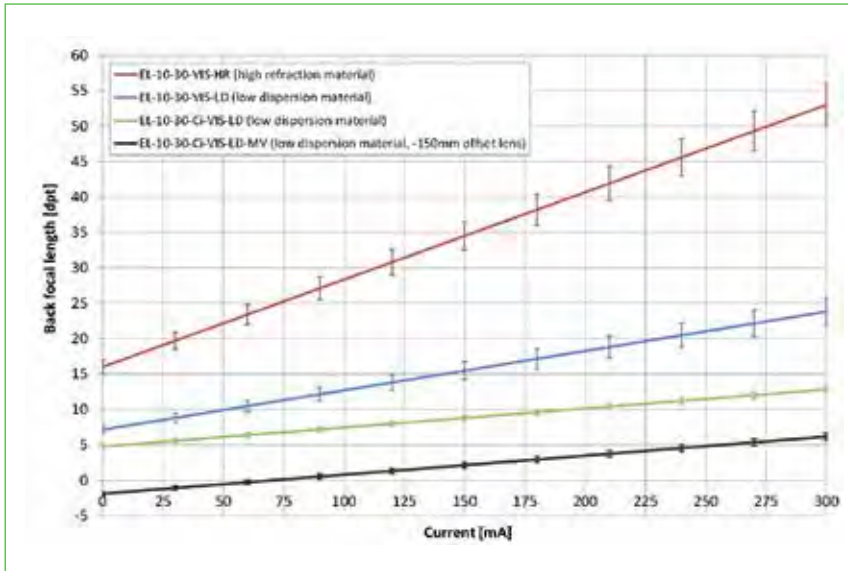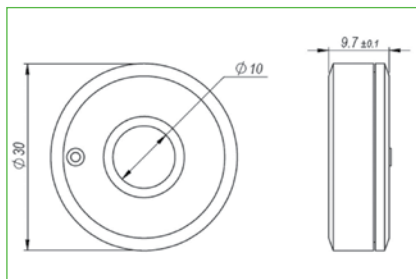
You can find more information and datasheets @ www.optotune.com

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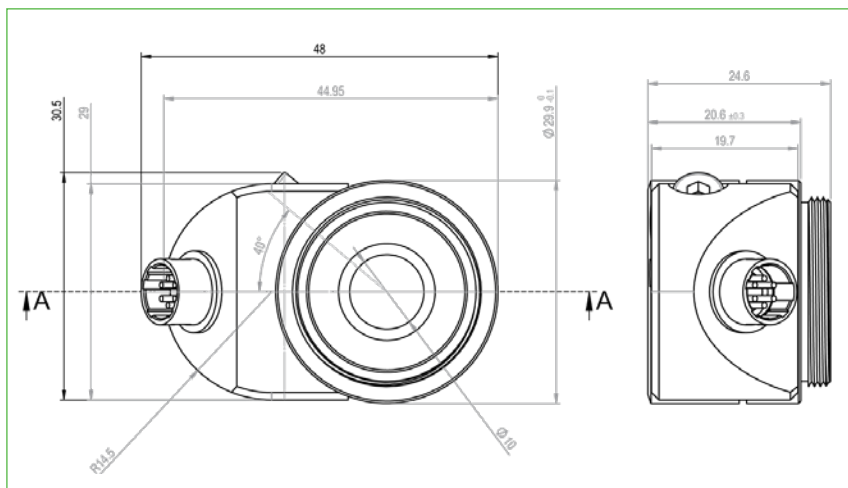Electrically Focus Variable Lens EL-16-40

By applying an electric current to this shape changing polymer lens, its optical power is controlled within milliseconds over a diopter range of -2 dpt to +3 dpt. A major hallmark of this shape changing lens is the clear aperture of 16mm. In order to achieve good repeatability and focus stability, Optotune integrated a temperature sensor allowing for in-situ compensation of temperature effects. The lenses are therefore ideally suited for a large variety of applications where imaging and a large beam diameter is of importance.

- Clear aperture: 16 mm
- Optical power: tuning range @ 30 °C with Optotune's Lens Driver 4: -2 to +3 dpt
- Repeatability: +/- 0.1 dpt
- Wavefront error (@525 nm, 0 mA) Optical axis vertical / horizontal: ~ 0.25 / ~0.5 λ RMS
- Lens type: plano-concave to plano-convex
- Cover glass coating: 420 to 950 nm
- Response time (typical at 30°C, 0 to +/- 250 mA step): 5 ms
- Settling time (typical at 30 °C, 0 to +/- 250 mA step): 25 ms
- Lifecycles (10% - 90% sinusoidal): >1.000.000.000
- Operating temperature: -20 to 65 °C
- Storage temperature: -40 to 85°C
- Nominal control current with Lens Driver 4: -250 to 250 mA
- Absolute max. control current: -500 to 500 mA
- Power consumption: 0 to 0.7 (nominal), 0 to 2.8 (absolute max.) W
- Motor coil resistance @ 30°C: 11 Ω
- Voltage for digital circuitry Vcc: 3.3 V

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Electrical Lens Driver

The Electrical Lens Driver is a simple-to-use tool for controlling Optotune Electrically Focus-Tunable Lenses or C-Mount Focus-Tunable Lenses. The USB powered Optotune Electrical Lens driver offers I2C sensor read-out for temperature compensation, in addition to rectangular, triangular, or sinusoidal drive frequencies from 0.2 –1000Hz. Additional on-board components include an 8-bit, 16MHz microcontroller with 32KB flash memory and a 6-way FPC connector. The driver is ideal for standalone use or for OEM integration.

- Standalone or for integration in OEM designs
- To control electrically tunable lenses of series EL-10-30
- Driver software download at www.optotune.com/downloads
- Control range of -200 to +200mA with 12-bit precision

Steuermodul Lens Driver 4 Electrical Lens Driver 4

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Electrical Lens Driver 4	G399579000
Electrical Lens Driver EL-E-4i	G399573000
Hirose cable CAB-6-100 (6 Pin, 100cm length)	G399574000
Hirose cable CAB-6-300 (6 Pin, 300cm length)	G399585000

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Profit from our 135 years of experience in the production of precision optical and mechanical components.

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*Optics production with CNC polishing
machine*

The LINOS Achromats and Lens Systems

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Our quality criteria:

- *High-quality materials from established manufacturers (Schott & Ohara)*
- *Extremely narrow focal length tolerances*
- *Minimal chromatic length aberration*
- *Minimal spherical aberration*
- *Minimal wavefront distortion (between $\lambda/4$ and $\lambda/10$)*
- *Compatible with LINOS and other bench systems*

Achromats & Lens Systems

Mounted or unmounted, coated or uncoated, with positive or negative focal lengths, LINOS achromatic lenses can fulfill every one of your optical requirements and exemplify our continued expertise in manufacturing complex optics.

LINOS achromatic lenses ensure consistently high imaging quality with resolutions that approach the theoretical diffraction limit. Laser centering and continuous, strict quality control guarantee the highest standards for surface form and alignment of the curvature center in every single achromatic lens that leaves our factory. Compare for yourself!

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Special Features:

All relevant structural data (radii, center thicknesses, materials) can be found in our WinLens optics software package. Broadband anti-reflective coatings and customer-specific coatings are available.

Ideal Areas:

Beam focusing, imaging systems, fiber collimation.

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

Content

Technical Information	286
Achromats	
Achromats VIS, Positive; Dia. 3 mm to 31.5 mm, Unmounted	288
Achromats VIS, Positive; Dia. 3 mm to 31.5 mm, Mounted	290
Achromats VIS, Positive; From 40 mm Diameter, Unmounted	292
Lens Systems	
Condenser Combination, $f=16$ mm, Crown Glass	293
Incident Light Microscope Objectives	293
HD Micro Objectives	294

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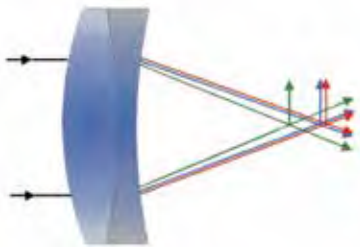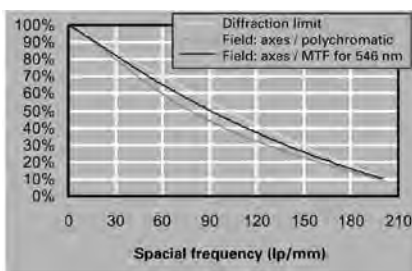
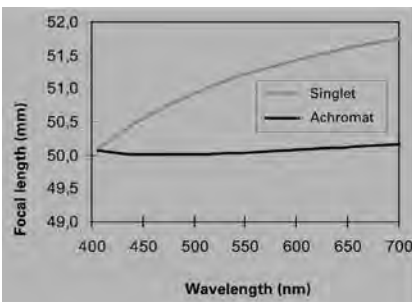
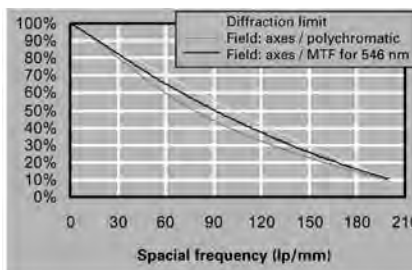
Comparison of the wavelength-dependent focal length
change of a single lens and an achromat

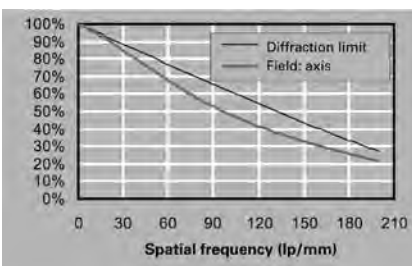Poly- and monochromatic MTF curve for achromat with
 $f = 200\text{ mm} / \varnothing 31.5\text{ mm}$

Polychromatic MTF curve for achromat with
 $f=120\text{ mm} / \varnothing 25.4\text{ mm}$

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An achromat with a positive refractive power consists of a convex crown glass lens with low dispersion (color splitting) and a concave flint glass lens with a high dispersion.

The lenses are matched such that the variation of the focus position of the overall system is minimized for two selected wavelengths. A modern chromatic (color) correction of the focus position is calculated for the Fraunhofer lines C' (479.9914 nm) and F' (643.8469 nm). The remaining variation of the focus position (secondary spectrum), e.g. to the e-line (546.074 nm), is typically below 0.1% of the focal length.

In addition, the spherical aberration can be minimized by suitable geometry and material selections. LINOS achromats are especially well corrected against spherical aberration and are therefore very good for monochrome applications for optimum focusing and/or collimating with minimum wave front deformation, e.g. of laser light.

In the design of the LINOS achromats special attention was paid to the coma correction close to the axis (observance of the Abbe's sine conditions). This type of correction makes the system insensitive to slight tilting. This way, small, extended objects can be mapped very well.

Achromats are optimized for imaging remote objects. Therefore, they achieve their best imaging performance with light incidence parallel to the axis (e.g. focusing collimated laser radiation). Achromats are asymmetric; therefore one must observe a correct alignment in the beam path. Typically, the more strongly curved the surface faces are the more parallel the light is.

All LINOS achromats are laser-centered and therefore have an excellent alignment of the center of curvature (optical axis) to the edge of the achromat. In combination with high demands on the surface shape, this ensures a constant high imaging quality. This optical imaging performance is subject of strict quality control. Typical wave front deformations are in a range between $\lambda/4$ and $\lambda/10$.

The modulation transfer function (MTF) illustrates the high imaging quality. This type of graph shows the resolution capacity in dependence of the resolution of the object or image structure (also see the chapter "Technical Information on Optics". The resolution is usually close to the theoretical diffraction limit.

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
LINOS achromats are standardly available with an ARB 2 VIS coating.

This coating is a hard and scratchproof anti-reflective coating.

Please refer to the reflection diagram for the characteristics of this coating.

One outstanding feature of this coating is the extrem high transmission of our achromats which is > 99 % for the stated wavelength range.

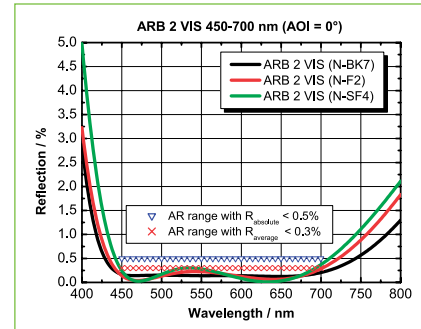Besides the ARB2 VIS for the visible spectral range, other coating as a ARB2 NIR (725-1050 nm) or a UV-VIS for a spectral wavelength range of 240-700 nm are available on request. More information about our coating competence are described in chapter „Thin Film Coatings“.

ARB2 Broadband-anti-reflective coating for VIS, Transmission for 450-700 nm > 99 %

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
C*	Optik-Ø (mm)	Sauberkeit ISO 10110	Ø Toleranz (mm)	dm Toleranz (mm)	Zentriergenauigkeit
C*	Ø optics (mm)	Surface Quality ISO 10110	Ø Tolerance (mm)	dm Tolerance (mm)	Centering Accuracy
1	3-6	3x0.1	-0.03	±0.2	better than 4'
2	8-12.5	3x0.16	-0.05	±0.2	better than 4'
3	18	5x0.16	-0.05	±0.2	better than 4'
4	22.4-25.4	3x0.16	-0.06	±0.2	better than 4'
5	31.5	5x0.25	-0.07	±0.2	better than 4'
6	40-80	5x0.4	-0.08	±0.2	better than 6'
7	100	5x0.63	-0.09	±0.2	better than 6'

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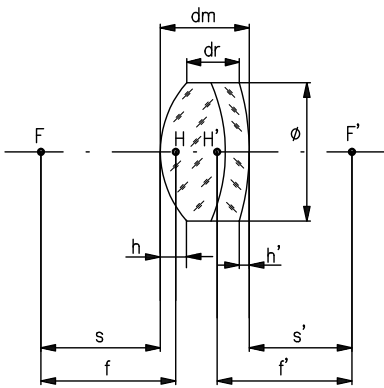Qioptiq uses glass from established manufacturers in the production of precision optics components. This ensures the consistent high quality of our lenses, achromatic lenses and plano-optics.

f focal length
s object distance
s' image distance
Ø lens diameter
dm center thickness
dr edge thickness
h primary vertex (h=dm-dr-h')
h' sagittal height secondary vertex
F, F' focal points
H, H' principal points

Tolerances:
Focal length f: ±1%
Image distance s': ±1%

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Achromats VIS, Positive; Dia. 3 mm to 31.5 mm, Unmounted

- Minimal longitudinal chromatic aberration
- Minimal spherical aberration
- Minimal wavefront distortion
- Tight focal length tolerance
- Broadband anti-reflection coated
- Cemented doublets
- Unmounted
- Broadband anti-reflection coating ARB2 VIS
- Damage thresholds > 2 kW/cm² cw (488/514 nm)
- Damage thresholds > 200 mJ/cm² for pulses of 10 ns (1064 nm)

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
All relevant specifications of the lenses, such as materials, radii and center thicknesses, are included in the database of our WinLens optical design software. A freeware version of this program is available for download at winlens.de.

Achromate VIS, Positiv, Durchmesser 3 - 31.5 mm, ungefasst Achromats VIS, Positive; Dia. 3 mm to 31.5 mm, Unmounted

Brennweite (mm)	Ø (mm)	dm (mm)	dr (mm)	h' (mm)	s' _{546 nm} (mm)	f' _{546 nm} (mm)	C ⁺	Part No.
Focal length (mm)	Ø (mm)	dm (mm)	dr (mm)	h' (mm)	s' _{546 nm} (mm)	f' _{546 nm} (mm)	C ⁺	Part No.
4	3	2.4	1.8	0.2	2.83	4	1	G32218000
5	3	3.1	2.6	0.2	3.6	5	1	G32219000
6	3	2.7	2.2	0.1	4.68	5.95	1	G322250000
8	6	5	3.7	0.4	5.48	8	1	G32220000
10	6	3	2.1	0.2	8.88	10.38	1	G322206000
13	6	4.9	4.1	0.2	10.53	13	1	G32222000
18	6	3.7	3.2	0.1	16.07	18	1	G32223000
16	8	3.5	2.4	0.2	14.53	16.32	2	G322207000
25	8	4.2	3.4	0.3	22.88	24.97	2	G322356322
30	8	4.2	3.5	0.2	27.86	29.95	2	G322357322
35	8	4.2	3.6	0.2	32.88	34.96	2	G322358322
40	8	4.2	3.7	0.15	37.88	39.96	2	G322359322
45	8	4.3	3.9	0.15	42.86	44.96	2	G322360322
60	8	4.4	4.1	0.1	57.8	59.94	2	G322362322
80	8	4.4	4.2	0.08	77.84	79.97	2	G322363322
100	8	4.5	4.3	0.06	97.73	99.9	2	G322364322

Brennweite (mm)	Ø (mm)	dm (mm)	dr (mm)	h' (mm)	s' _{546 nm} (mm)	f' _{546 nm} (mm)	C ⁺	Part No.
Focal length (mm)	Ø (mm)	dm (mm)	dr (mm)	h' (mm)	s' _{546 nm} (mm)	f' _{546 nm} (mm)	C ⁺	Part No.
150	8	4.5	4.4	0.04	147.75	149.92	2	G322366322
200	8	4.7	4.6	0.03	197.49	199.73	2	G322367322
300	8	4.7	4.6	0.02	297.61	299.86	2	G322368322
20	10	4.1	2.8	0.3	18.44	20.6	2	G322201000
25	12.5	5.5	4	0.3	22.38	25.33	2	G322284000
30	12.5	5	3.7	0.3	27.59	30.24	2	G322285000
35	12.5	5	3.9	0.2	32.39	35.04	2	G322286000
40	12.5	6.3	5.1	0.4	37.24	40.09	2	G322337000
45	12.5	5.5	4.4	0.3	42.51	45.02	2	G322338000
50	12.5	5	4.1	0.3	47.65	50.01	2	G322339000
60	12.5	5.3	4.5	0.3	57.69	59.99	2	G322341000
80	12.5	5.5	5.1	0.2	77.57	79.94	2	G322343000
100	12.5	5.3	4.8	0.2	97.81	99.99	2	G322345000
120	12.5	5.8	5.4	0.12	117.93	119.91	2	G322320322
150	12.5	5.8	5.5	0.1	147.42	149.9	2	G322325322
200	12.5	5.5	5.3	0.08	197.55	199.87	2	G322327322
300	12.5	4	3.8	0.1	297.57	299.85	2	G322348000
40	18	7	4.5	0.6	36.04	40.16	3	G322209000
50	18	6.5	4.6	0.3	45.95	50.04	3	G322265000
60	18	6	4.4	0.2	56.33	60.11	3	G322266000
80	18	5.8	4.6	0.2	76.62	80.13	3	G322210000
100	18	5.8	4.8	0.1	96.43	100.11	3	G322236000
150	18	6.5	5.9	0.18	146.94	149.83	3	G322331322
200	18	6.5	6.1	0.15	197.07	199.82	3	G322332322
50	22.4	7.5	5.2	0.4	45.76	50.04	4	G322321000
60	22.4	7	5.1	0.4	56.74	60.47	4	G322322000
80	22.4	7	5.1	0.6	77.05	80.34	4	G322323000
100	22.4	9	7.5	0.5	95.89	100.16	4	G322324000
140	22.4	8	6.9	0.4	136.2	140.07	4	G322326000
160	22.4	7	6.5	0.3	156.6	159.8	4	G322333322
200	22.4	6	5.2	0.2	197.38	200.26	4	G322328000
400	22.4	7	6.6	0.12	396.5	399.6	4	G322334322
500	22.4	6	5.7	0.1	495.99	499.43	4	G322329000
60	25.4	8.9	6.2	0.7	55.47	60.1	4	G322306322
80	25.4	7	4.6	0.8	77.05	80.34	4	G322307322
100	25.4	9	7.1	0.6	95.89	100.16	4	G322308322
120	25.4	7	5.4	0.5	116.77	120.1	4	G322309322
140	25.4	8	6.6	0.5	136.2	140.07	4	G322351000
160	25.4	8.6	7.4	0.4	155.95	160	4	G322352000
200	25.4	6	5	0.3	197.38	200.26	4	G322353000
300	25.4	7	6.4	0.2	296.5	299.66	4	G322336322
400	25.4	7	6.5	0.15	396.44	399.61	4	G322340322
500	25.4	6	5.5	0.1	495.99	499.43	4	G322355000
80	31.5	10.5	6.7	1.2	74.84	80.1	5	G322287322
100	31.5	9.5	6.5	1	95.34	100.09	5	G322288322
120	31.5	9.5	7	0.8	115.6	119.91	5	G322269322
140	31.5	8	5.9	0.7	136.2	140.07	5	G322239322
150	31.5	8.6	6.6	0.7	145.85	150.03	5	G32227322
160	31.5	8.6	6.7	0.6	155.95	160	5	G322270322
180	31.5	9	7.1	0.63	175.84	180.1	5	G322342322
200	31.5	7	5.5	0.4	197.21	200.41	5	G322271322
250	31.5	7	5.8	0.4	247.42	250.45	5	G322272322
300	31.5	7	6	0.3	297.19	300.23	5	G322273322
400	31.5	7	6.3	0.2	397.05	400.08	5	G322275322

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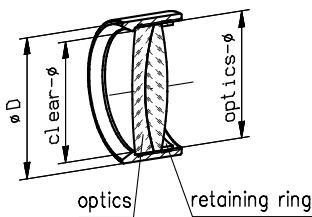
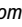
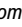
Achromats VIS, Positive; Dia. 3 mm to 31.5 mm, Mounted

- Minimal longitudinal chromatic aberration
- Minimal spherical aberration
- Minimal wavefront distortion
- Tight focal length tolerance
- Broadband anti-reflection coated with ARB2 VIS
- Achromats in Nanobench  and Microbench  mounts fitting directly to these LINOS systems

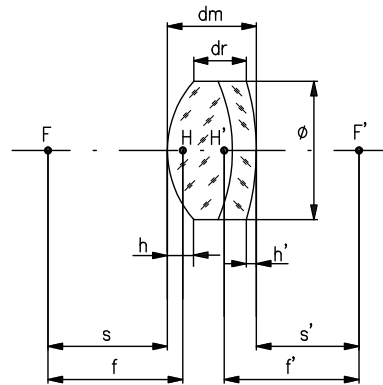- Cemented doublets
- Damage thresholds > 2 kW/cm² cw (488/514 nm)
- Damage thresholds > 200 mJ/cm² for pulses of 10 ns (1064 nm)

f	focal length
s	object distance
s'	image distance
\varnothing	lens diameter
dm	center thickness
dr	edge thickness
h	primary vertex ($h=dm-dr-h'$)
h'	sagittal height secondary vertex
F, F'	focal points
H, H'	principal points

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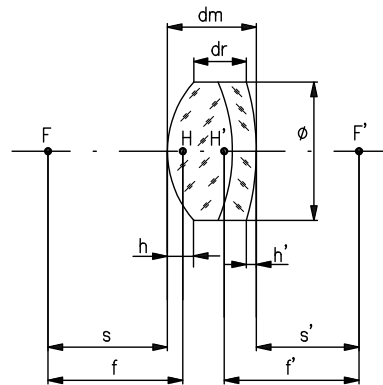
Achromats VIS, Positive; From 40 mm Diameter, Unmounted

- Minimal longitudinal chromatic aberration
- Minimal spherical aberration
- Minimal wavefront distortion
- Tight focal length tolerance
- Broadband anti-reflection coated with ARB2 VIS
- Cemented doublets
- Damage thresholds > 2 kW/cm² cw (488/514 nm)
- Damage thresholds > 200 mJ/cm² for pulses of 10 ns (1064 nm)


f focal length
s object distance
s' image distance
Ø lens diameter
dm center thickness
dr edge thickness
h primary vertex (h=dm-dr-h')
h' sagittal height secondary vertex
F, F' focal points
H, H' principal points

Tolerances:
Focal length f: ±1%
Image distance s': ±1%

Achromate VIS, Positiv, ab 40 mm Durchmesser, ungefasst Achromats VIS, Positive From 40 mm Diameter, Unmounted

Brennweite (mm)	Ø (mm)	dm (mm)	dr (mm)	h' (mm)	s' _{546 nm} (mm)	f' _{546 nm} (mm)	f' _{1064 nm} (mm)	⊙*	Part No.
Focal length (mm)	Ø (mm)	dm (mm)	dr (mm)	h' (mm)	s' _{546 nm} (mm)	f' _{546 nm} (mm)	f' _{1064 nm} (mm)	⊙*	Part No.
90	40	14.3	9.8	0.8	81.78	89.98	90.83	6	G322389000
120	40	13	9.7	0.9	113.6	119.99	120.96	6	G322388000
160	40	11.7	9	0.8	154.64	160.21	161.35	6	G322384000
180	40	10.8	8.2	0.4	174.92	181.66	183.18	6	G322246000
200	40	10	7.7	0.4	195.34	201.26	203.02	6	G322293000
250	40	9.5	7.6	0.6	245.52	249.91	251.7	6	G322294000
300	40	11	9.4	0.5	294.78	300.09	302.77	6	G322295000
80	50	20	11.9	1.7	69.05	80.14	80.65	6	G322301000
100	50	17.5	11.2	1.4	92.03	101.57	102.06	6	G322302000
120	50	15.5	10.3	1.1	112.88	121.36	121.82	6	G322303000
200	50	12.5	8.9	0.6	193.7	201.27	203.05	6	G322304000
300	50	12	9.6	0.7	294.31	300.18	302.86	6	G322305000
160	50.8	13.5	8.6	1.5	152.76	159.49	160.72	6	G322310000
250	50.8	11.5	8.4	1	244.65	250.1	251.9	6	G322311000
400	50.8	10.5	8.6	0.6	395.59	400.33	403.14	6	G322312000
150	63	21	13.9	1.3	138.3	150.44	151.51	6	G322383000
500	63	13	10.7	0.7	494.65	500.39	503.9	6	G322230000
800	63	13	11.5	0.5	796.93	802.79	808.31	6	G322232000
160	80	29	18.8	2.1	144.45	160.1	160.89	6	G322267000
310	80	18.5	12.3	1.9	301.5	310.35	312.69	6	G322278000
500	80	16	12.2	1.2	493.2	500.64	504.17	6	G322279000
500	100	22	16	1.9	488.87	499.43	503.04	7	G322316000
1000	100	21.5	18.5	0.6	988.52	1000.36	1007.49	7	G322313000

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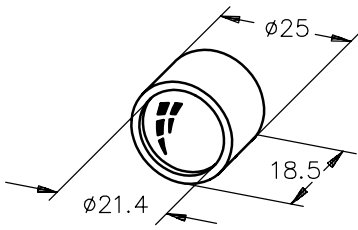
Condenser Combination, f=16 mm, Crown Glass

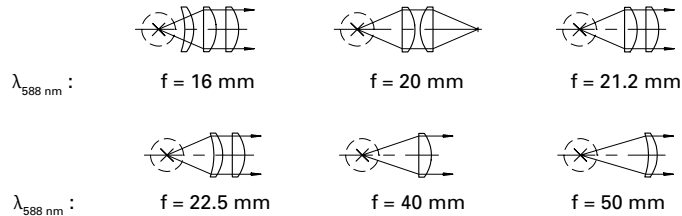- For uniform illumination
- For use as collimators in large focal ratio illumination systems
- Supplied in CL-mounts, Ø 25 mm, 21.4 mm clear-aperture
- Microbench-compatible
- Re-arrangeable lens elements (see illustrations; all tools required are included)
- Consisting of: 2 plano-convex lenses f=40 mm (G31 2314 000), 1 concave-convex lens f=50 mm (G31 5505 000)
- For customized optics and higher volumes please contact our sales team for your individual quotation.

Kombinationskondensor, f = 16 mm aus Kronglas Condenser combination, f = 16 mm, crown glass

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Condenser combination, f = 16 mm, crown glass	G063010000

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Incident Light Microscope Objectives

- For use with incident illumination; without cover glass correction
- Also suitable for laser beam expanding
- Achromatic correction
- Parfocal length 45 mm
- RMS thread (W0.8"×1/36")
- Thread length 4.5 mm

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HD Micro Objectives

Owed to many years of experience in the field of Chip-On-The-Tip technology, Qioptiq was able to advance the development and production of new objectives for small image sensors.

Significant features of the 1/10"-FALCON series are the high MTF and the unvignetted image.

All types have an integrated infra-red-absorbing filter and a lithographically generated aperture.

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Our quality criteria:

- *High-quality glass materials from established manufacturers*
- *Many years of experience in precision optics production*
- *The narrowest possible production tolerances*

LINOS Plano Optics

Qioptiq is known for high-quality plane plates from optical glass, fused silica and sapphire, as well as a variety of prisms, filters, beamsplitters, reticles and dispersion plates. We can offer support for all your applications that require precise planarity and flawless surfaces.

High-quality materials ensure the best possible transmission and consistent, reliable quality standards in all Qioptiq products.

In production, we place a high value on the narrowest possible angular tolerances and the best parallelism, integrity and fit. Qioptiq plano optics and their coatings are manufactured using the latest machines and are subject to constant, meticulous quality control. Qioptiq stands for lasting precision!

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Special Features:

Many components are available in mounts compatible with LINOS bench systems. Prisms can also be found in the Prism Wizard of the WinLens 3D program.

Ideal Areas:

Beam deflection, Beam splitting, Measurement tasks, View panels, and more.

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Content

<i>Reflective Curves of Coated Plane Plates</i>	298
<i>Price List for Standard Coatings</i>	299
Plane Plates	
<i>N-BK7 Plane Plates</i>	300
<i>UV-grade Fused Silica Plane Plates, Standard Quality</i>	301
<i>UV-grade Fused Silica Plane Plates, High Flatness</i>	302
<i>UV-grade Fused Silica Plane Plates, High Flatness, High Parallelism</i>	302
<i>Fused Silica Brewster Plates</i>	303
<i>Shear Plates</i>	303
Prisms	
<i>90°-Prisms (isosceles)</i>	304
<i>90°-Prisms (Silver- and Anti-Reflection Coated)</i>	305
<i>90°-Porro Prisms (Isosceles)</i>	306
<i>Penta Prisms</i>	306
<i>60°-Prisms (Equilateral)</i>	307
Filters	
<i>UV Colored Glass Filters</i>	308
<i>Blue Colored Glass Filters</i>	309
<i>Green Colored Glass Filters</i>	310
<i>Colored Glass Filters, Different Edge-Positions</i>	311
<i>Color Filters DT</i>	312
<i>Color Filters DC</i>	314
<i>Neutral-Density Filters</i>	314
<i>Neutral Wedges</i>	316
<i>Heat-Absorbing Filters</i>	316
<i>Heat-Protection Filters CALFLEX™ B1/K1</i>	317
<i>Heat-Protection Filters CALFLEX™ 3000 SP</i>	317
<i>Heat-Protection Filters CALFLEX™ C</i>	318
<i>Cold Light Mirror KS 93 / 45°</i>	318
<i>NIR-Blocking Filters CALFLEX™ X</i>	319
<i>UV-Blocking Filters UV-B</i>	319
Beamsplitter	
<i>Beamsplitter Cubes UV</i>	320
<i>Beamsplitter Cubes VIS</i>	321
<i>Beamsplitter Cubes NIR</i>	322
<i>Beamsplitter Plates UV</i>	323
<i>Beamsplitter Plates VIS-NIR</i>	324
<i>Beamsplitter Plates NIR</i>	326
<i>Beamsplitter Plates IR</i>	327
<i>Beamsplitter Plates with Different Splitting Ratios</i>	328
Reticles and Dispersion Plates	
<i>Reticles</i>	329
<i>Surface Dispersion Plates (Ground Glass)</i>	330

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Reflective Curves of Coated Plane Plates

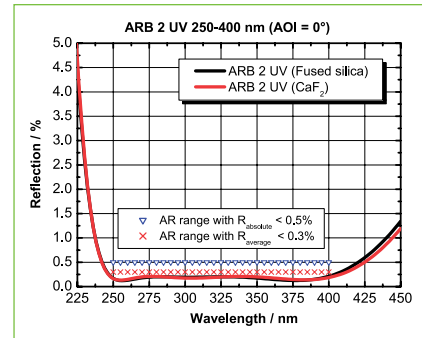
All crown and fused silica plane plates of the following pages are also available as coated standard items with rapid delivery.

UV Broadband Anti-Reflective Coating ARB2 UV

- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the UV spectral range*
- *Guaranteed residual reflection $< 0.5\%$ for angles of incidence $0^\circ \leq \text{AOI} \leq 15^\circ$ in the AR range*
- *Typical average residual reflection $< 0.3\%$ for angles of incidence $0^\circ \leq \text{AOI} \leq 15^\circ$ in the AR range*
- *Damage threshold $H_{\infty} > 2 \text{ J/cm}^2$ at 266 nm with 9 ns laser pulses (s-on-1) at 10 Hz*
- *AR range, typical: 250-400 nm*
- *Design adaptable in the spectral range of $245 \text{ nm} \leq \lambda \leq 450 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, CaF_2*

Ordering Information:

To order plane plates with ARB2 UV on both sides simply change the last three numbers of the order number to 522.

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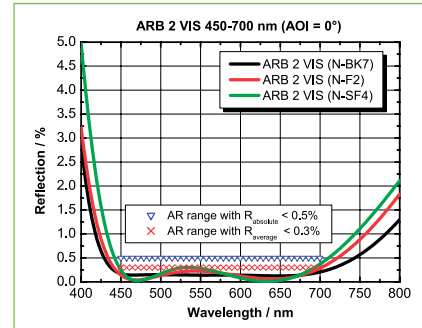
Broadband Anti-Reflective Coating ARB 2 VIS

- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the visible spectral range*
- *Guaranteed residual reflection $< 0.5\%$ for angles of incidence $0^\circ \leq \text{AOI} \leq 15^\circ$ in the AR range*
- *Typical average residual reflection $< 0.3\%$ for angles of incidence $0^\circ \leq \text{AOI} \leq 15^\circ$ in the AR range*
- *Damage threshold $H_{\infty} > 10 \text{ J/cm}^2$ at 532 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *AR range typical: 450-700 nm*
- *Design adaptable in the spectral range of $350 \text{ nm} \leq \lambda \leq 950 \text{ nm}$ and for higher angles of incidence*
- *Substrates: N-BK7, fused silica or similar*

Ordering Information:

Simply change the last three numbers of the order number to 322.

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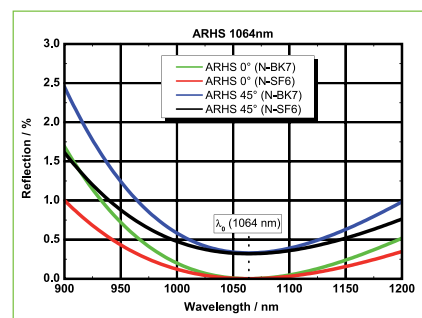
High-power narrowband antireflective coating ARHS

- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for 1064 nm*

Ordering Information:

Simply change the last three numbers of the order number to 329.

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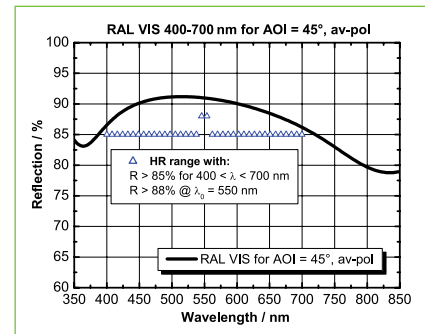
Reflective aluminum coating RAL VIS

- *Standard aluminium reflective coating for mirrors in the VIS and IR spectral range up to 3 µm*
- *High reflection within a broad spectral and angle-of-incidence range*
- *Guaranteed reflection > 85 % in the range of 400 nm ≤ λ ≤ 700 nm, > 88 % at λ₀ = 550 nm for AOI = 45°, unpolarized*
- *Damage threshold H_∞ > 1.5 J / cm² at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *Design adaptable in the spectral range of 350 nm ≤ λ ≤ 3000 nm and for higher angles of incidence*
- *Substrates: optical glasses, ceramic glasses, Zerodur, Borofloat, Pyrex, or similar*

Ordering Information:

Simply change the last three numbers of the order number to 331.

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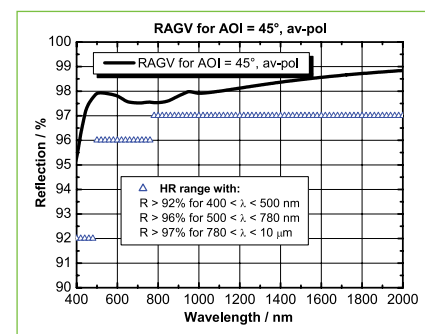
Front surface silver coating RAGV

- *Front surface mirror silver coating for the VIS and IR spectral range up to 12 µm*
- *Constant high reflection within a broad spectral and angle-of-incidence range without absorption bands*
- *Guaranteed reflection (0° ≤ AOI ≤ 60°, unpolarized) > 96 % in the range of λ > 500 nm*
- *Damage threshold H_∞ > 1.5 J/cm² at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *Substrates: optical glasses, ceramic glasses, Zerodur, Borofloat, Pyrex, or similar*

Ordering Information:

Simply change the last three numbers of the order number to 400.

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Note:

If fused silica is used as a substrate material, the curves of antireflection coatings designed for N-BK7 are slightly modified.

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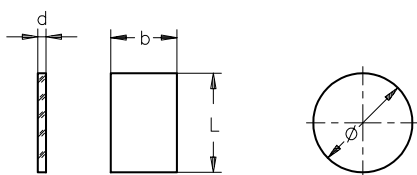
N-BK7 Plane Plates

- Ideal substrates for planar optical components, such as plane mirrors, filters or beamsplitter plates
- Ideal for use as windows in optical systems
- Surfaces uncoated

Rectangular type and circular type

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For higher quantities, our plane plates are also available with the following standard, economically priced anti-reflective coatings:

- ARB 2 VIS (450-700 nm), order code -322
- ARB 2 NIR (725-1050 nm), order code -525 (on request)
- ARHS 1064 (Nd:YAG center wavelength at 1064 nm), order code -329

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When placing an order or requesting a quote, simply replace the "000" in the order number with the order code for the desired coating. For more information please look at the beginning of this chapter respectively chapter "Thin Film Coatings."

Planplatten aus N-BK7 Plane plates fabricated from N-BK7

Größe bxL oder Ø (mm)	Dicke d (mm)	Toleranz d (mm)	Toleranz b (mm)	Toleranz L (mm)	Planität beidseitig (λ)	Toleranz Ø (mm)	Keilwinkel γ	Part No.
Dimensions bxL or Ø (mm)	Thickness d (mm)	Tolerance d (mm)	Tolerance b (mm)	Tolerance L (mm)	Flatness Both Surfaces (λ)	Tolerance Ø (mm)	Wedge Angle γ	Part No.
15x20	2	±0.2	-0.11	-0.13	1		2'	G390136000
20x30	2.5	±0.3	-0.13	-0.13	2		2'	G390137000
30x50	4	±0.2	-0.13	-0.16	2		3'	G390138000
Ø 10	2	±0.1			1/4	-0.09	2'	G390130000
Ø 12.7	2	±0.1			1/4	-0.11	1'	G390068000
Ø 16	2	±0.1			1/4	-0.11	2'	G390131000
Ø 22.4	3	±0.1			1/4	-0.13	3'	G390132000
Ø 25.4	5	±0.1			1/4	-0.15	1'	G390069000
Ø 31.5	3	±0.1			1/2	-0.16	3'	G390043000
Ø 40	6	±0.1			1/2	-0.16	3'	G390133000
Ø 50	6	±0.1			1/2	-0.16	3'	G390134000
Ø 63	6	±0.1			1	-0.19	3'	G390135000
Ø 80	10	±0.2			1	-0.19	3'	G390020000
Ø 100	10	±0.15			2	-0.22	3'	G390044000
Ø 150	15	±0.2			3	-0.25	5'	G390110000
Elliptical 22.4x31.5	3.5	±0.2			1/2		3'	G390139000

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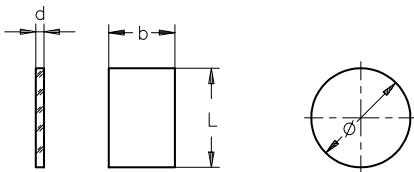
UV-grade Fused Silica Plane Plates, Standard Quality

- Ideal substrates for planar optical components, such as plane mirrors, filters, or beamsplitter plates
- Ideal for use as windows in optical systems
- Surfaces uncoated

Rectangular type and circular type

Planplatten aus Quarzglas mit Standardqualität Plane plates fabricated from UV-grade fused silica; standard quality

Größe bxL oder Ø (mm)	Dicke d (mm)	Toleranz d (mm)	Toleranz b (mm)	Toleranz L (mm)	Planität beidseitig (λ)	Toleranz Ø (mm)	Keilwinkel γ	Part No.
<i>Dimensions bxL or Ø (mm)</i>	<i>Thickness d (mm)</i>	<i>Tolerance d (mm)</i>	<i>Tolerance b (mm)</i>	<i>Tolerance L (mm)</i>	<i>Flatness Both Surfaces (λ)</i>	<i>Tolerance Ø (mm)</i>	<i>Wedge Angle γ</i>	<i>Part No.</i>
15x20	2	±0.2	-0.11	-0.13	1		3'	G390198000
20x30	2.5	±0.2	-0.13	-0.13	2		3'	G390158000
30x50	4	±0.2	-0.13	-0.16	3		3'	G390199000
Ø 10	2	±0.1			1/4	-0.09	3'	G390150000
Ø 12.7	2	±0.1			1/4	-0.11	1'	G390166000
Ø 16	2	±0.1			1/4	-0.11	3'	G390192000
Ø 22.4	3	±0.2			1/4	-0.13	3'	G390151000
Ø 25.4	5	±0.1			1/4	-0.15	1'	G390167000
Ø 31.5	3	±0.2			1/2	-0.16	3'	G390152000
Ø 40	4	±0.2			1/2	-0.16	3'	G390153000
Ø 50	5	±0.2			1/2	-0.16	3'	G390154000
Ø 80	8	±0.2			1	-0.22	3'	G390156000
Ø 100	10	±0.2			2	-0.25	3'	G390157000
Elliptical 22.4x31.5	3.5	±0.2			1/2		3'	G390159000

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The mounts you need for mounting plano optics in our bench systems are listed in chapter "Mounts and Posts", "Microbench", and "Tube System":

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UV-grade Fused Silica Plane Plates, High Flatness

- Ideal substrates for planar optical components, such as plane mirrors, filters, or beamsplitter plates
- Ideal for use as windows in optical systems
- Superior quality plates with high flatness of $\lambda/10$ on one surface (marked)
- 4 diamond polishing: better than 1 nm rms
- Surfaces uncoated

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UV-grade Fused Silica Plane Plates, High Flatness, High Parallelism

- Ideal substrates for planar optical components, such as plane mirrors, filters, or beamsplitter plates
- Ideal for use as windows in optical systems
- Superior quality circular plates with high flatness and high parallelism
- Surfaces uncoated

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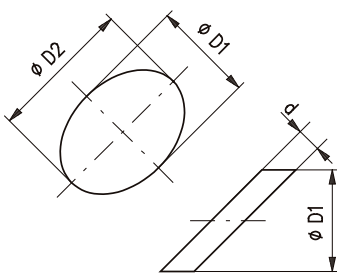
For plane mirrors in higher quantities, we offer the following standard, economically priced metallic mirror coatings for our plane plates:

- Aluminum mirror coating: RAL (400-3000 nm), order code -331
- Silver mirror coating: RAGV (450-3000 nm), order code -400

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For more information please look at the beginning of this chapter.

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
Fused Silica Brewster Plates

- To polarize light partially
- Ideal low-loss windows for use with polarized light
- Elliptically cut, bevelled edges
- Surfaces uncoated
- For customized optics and higher volumes please contact our sales team for your individual quotation.

Brewster-Platten aus Quarzglas

Brewster-plates fabricated from fused silica

Größe Ø D1 x Ø D2 (mm)	Dicke d (mm)	Toleranz d (mm)	Planität beidseitig (λ)	Toleranz Ø (mm)	Parallelität	Sauberkeit nach ISO 10110	Part No.
Dimensions Ø D1 x Ø D2 (mm)	Thickness d (mm)	Tolerance d (mm)	Flatness Both Surfaces (λ)	Tolerance Ø (mm)	Parallelism	Surface Quality	Part No.
12.7 x 22.5	4	± 0.2	1/10	-0.11	3'	3x0.063	G390215000
15.0 x 26.5	5	± 0.2	1/10	-0.11	3'	5x0.063	G390212000
25.4 x 45.0	6	± 0.2	1/10	-0.13	3'	5x0.063	G390211000

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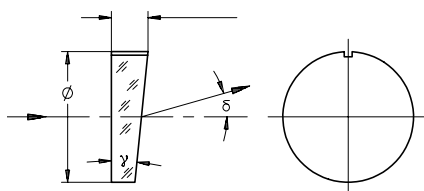
Shear Plates

- Wedge plate with high accuracy
- For collimation testing of laser beams
- For producing minimum angular or linear beam displacements
- Useful aid for equalizing minimum optical path differences
- N-BK7
- Surfaces uncoated
- λ/4 flatness on both faces

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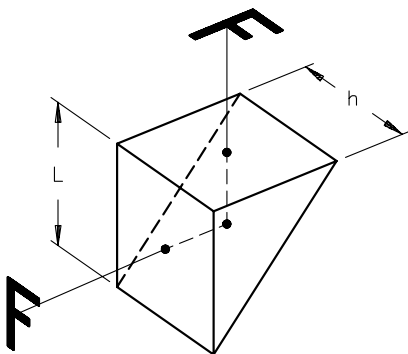90°-Prisms (isosceles)

- Deflect light rays
- Easy alignment
- Uncoated
- Please look at chapter Thin Film Coatings regarding customized coatings.
- Standard coating ARB2 for the visible spectral range or RAL/RAGV coatings are available for higher quantities on request.

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Making it easy to simulate and optimize off-axis systems: The Prism Wizard in our WinLens 3D optical design software is a user-friendly tool for implementation and modification of our standard prisms in models of optical systems. Try it! A freeware version of the program is available for download at www.winlens.de.

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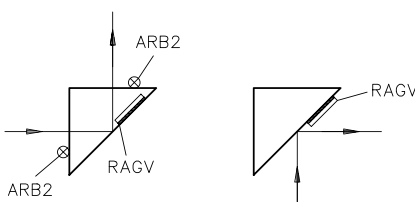
90°-Prisms (Silver- and Anti-Reflection Coated)


- Deflect light rays
- Easy to align
- To be used as deflection mirror or reflection prism
- Higher acceptance angle than uncoated prisms
- Hypotenuse RAGV (silver) coated ($R > 96\%$ for $\lambda = 500 - 1200\text{ nm}$)
- Angular deflection $\pm 10\text{ arcmin}$
- Prism tolerances max. 8 arcmin
- Made of N-BK7

To be used as deflection mirror or reflection prism

90°-Prismen (silberbeschichtet u. antireflexbeschichtet) 90°-Prisms (silver- and anti-reflection coated)

Abmessungen (mm)	Toleranz h (mm)	Toleranz L (mm)	Planität Hypotenuse (λ)	Planität Kathete (λ)	Part No.
Dimensions (mm)	Tolerance h (mm)	Tolerance L (mm)	Flatness Hypotenuse (λ)	Flatness Short Faces (λ)	Part No.
10x10	-0.09	-0.15	1/10	1/2	G339952000
12.5x12.5	-0.11	-0.15	1/10	1/2	G339959000
20x20	-0.13	-0.21	1/4	1	G339953000
25x25	-0.13	-0.21	1/4	1	G339955000
30x30	-0.13	-0.21	1/4	1	G339956000

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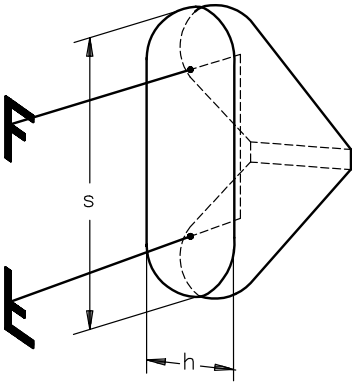
90°-Porro Prisms (Isosceles)

- 180° beam deflection, image-inverting
- Less alignment complexity than necessary with mirrors
- Porro principle
- N-BaK4 glass
- Uncoated

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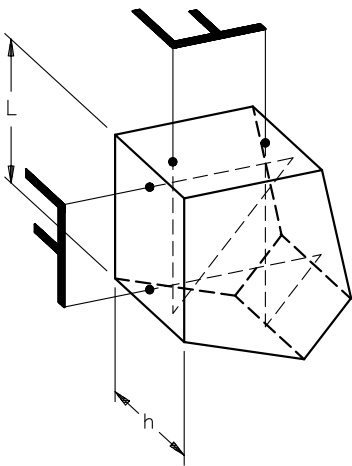
Penta Prisms

- Deflect beams through 90°
- No image reversal
- Beam deflection with different accuracy
- N-BK7 glass
- Coated reflective surfaces (RAGV)
- Entrance and exit faces broadband AR-coated (ARB 1)

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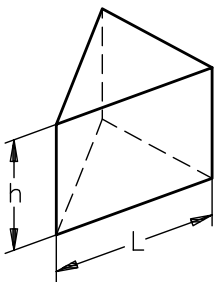60°-Prisms (Equilateral)

- Spectrally disperse light
- Fused silica versions for UV applications available
- Angular dispersion $\Delta\epsilon' = \epsilon'_{F'} - \epsilon'_{C'}$ ($\epsilon'_{F'}$ is the exit angle for $\lambda = 480$ nm, $\epsilon'_{C'}$ is the exit angle for $\lambda = 643.8$ nm)
- Uncoated

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Compatible mounts for prisms and beamsplitters are available in the chapters "Microbench", "Nanobench", "Tube Mounting System C", "Mounts and Posts" and "Positioning Systems".

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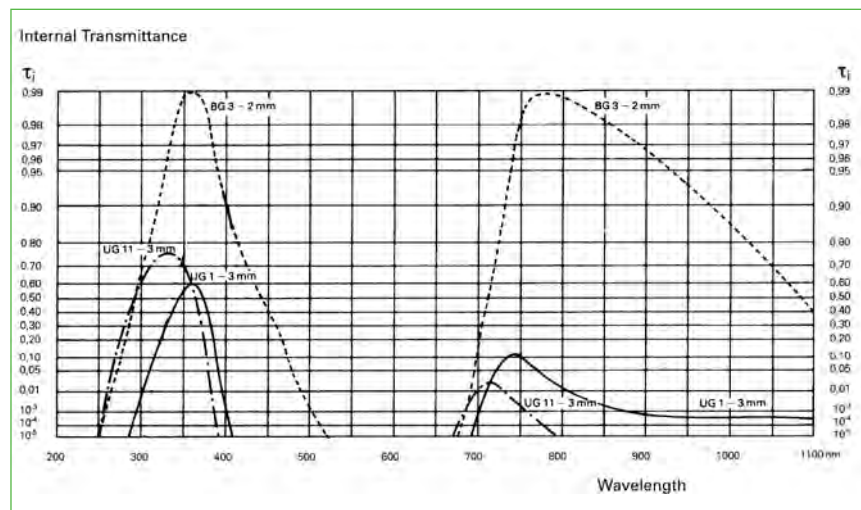
☒ In mount fitting directly to Microbench, see chapter "Mounts, Posts and Holders", section "Lens Mounts" for mount dimensions.

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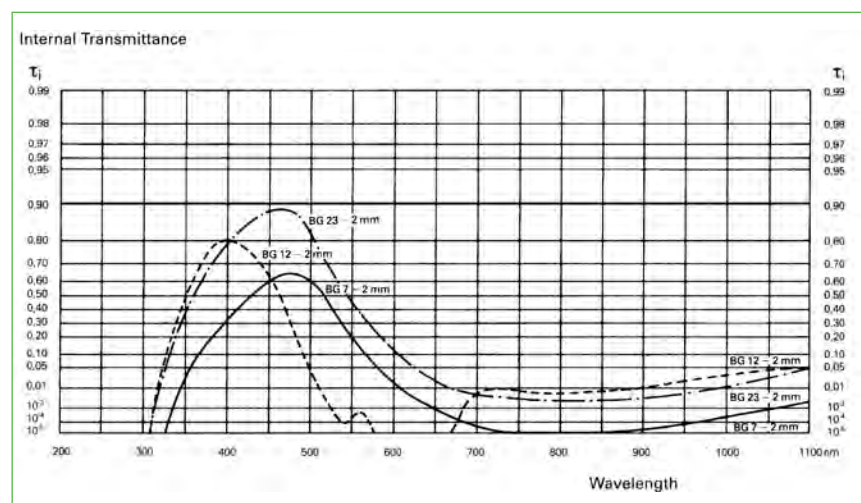
Blue Colored Glass Filters

- Volume filters fabricated from colored glass
- Mounted and unmounted versions available
- Mounted types fit to LINOS Microbench ☒

Tolerances:

- Ø 22.2 mm: ±0.2 mm; Ø 50 mm: +0/-0.6 mm
- Thickness: ±0.2 mm
- Bubbles: 1/5x0.25
- Striations: 2/03
- Surface Finish: 5/3x0.4
- Surface Parallelism: 15'

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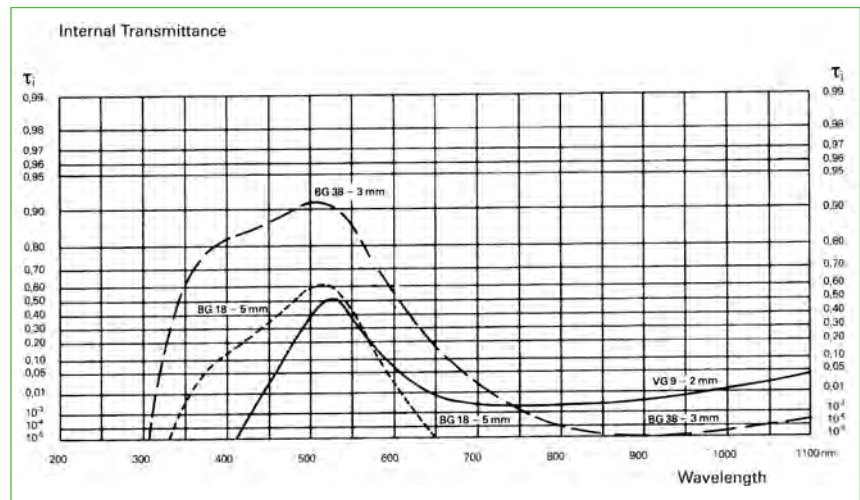
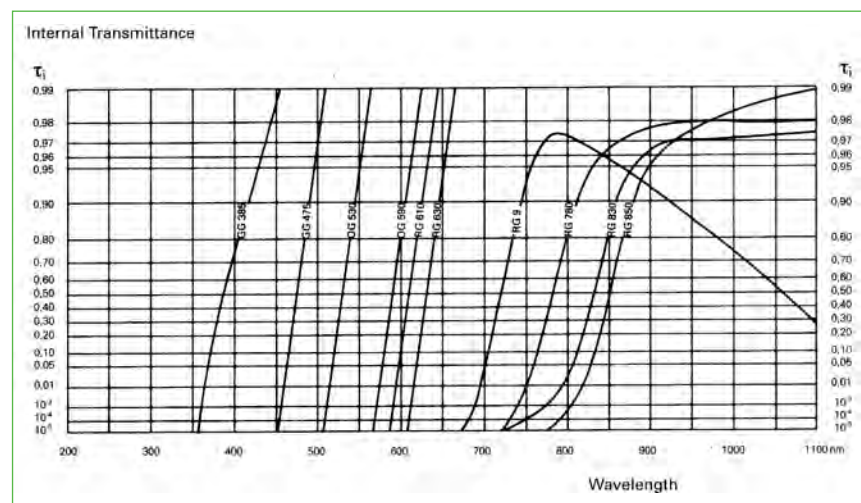
Colored Glass Filters, Different Edge-Positions

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- Mounted and unmounted versions available
- Mounted types fit to LINOS Microbench ☒

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- Tolerances:
- Ø 22.2mm: ± 0.2 mm; Ø 50 mm: +0/-0.6
 - Thickness: ±0.2 mm
 - Bubbles: 1/5x0.25
 - Striations: 2/03
 - Surface Finish: 5/3x0.4
 - Surface Parallelism: 15'

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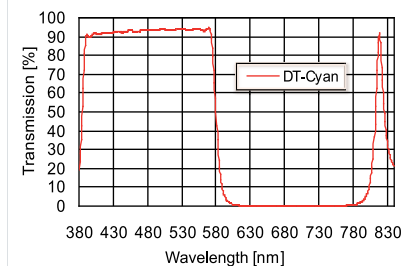
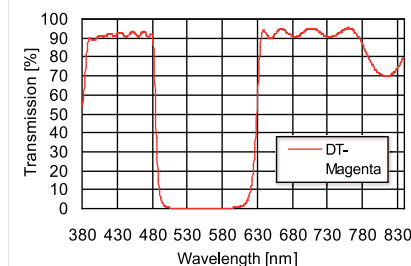
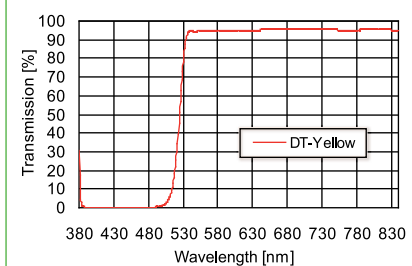
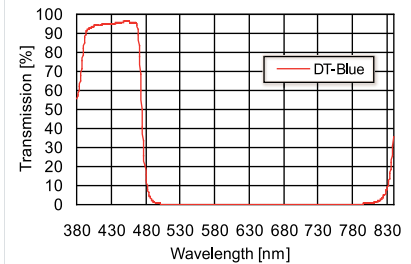
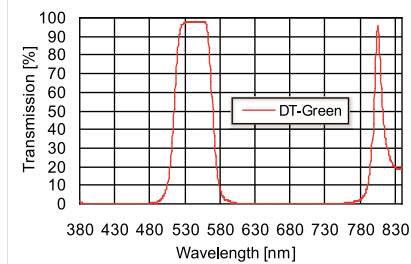
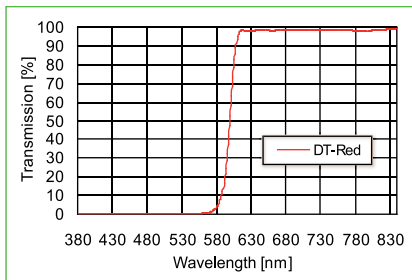
Color Filters DT

- Absorption free interference systems
 - Reflectance / transmittance of selected parts of VIS with high efficiency
 - Edge filters made to small tolerances for angles of incidence of 0°
 - Extremely high color saturation
 - Hard, dense dichroic coatings
 - Manufactured by patented "Sputter" technique
 - High temperature stability of spectral features
 - Heat-resistant up to 400°C
 - No aging or fading
 - High mechanical resistance
- Thickness of all filters: 1.1 mm
 - Special sizes up to 160 mm x 110 mm on request.

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
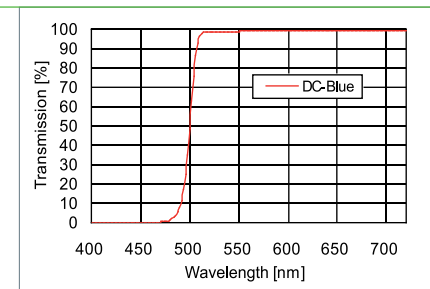
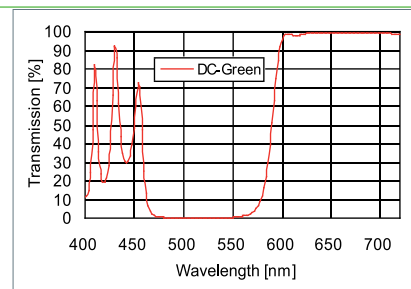
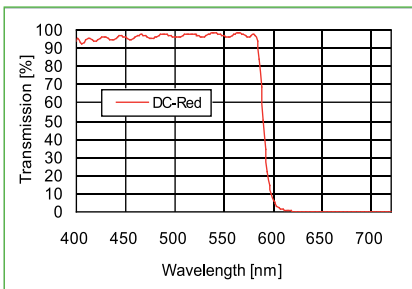
Color Filters DC

- Absorption free interference systems
- Nearly no-loss, diversion or mixture of basic colors blue, green and red
- For angles of incidence of 45°
- Heat-resistant up to approx. 300 °C
- Thickness of all filters: 1.1 mm
- Special sizes up to 160 mm x 110 mm on request.

Farbfilter DC Color Filters DC

Typ	Abmessungen (mm)	Part No.
Type	Dimensions (mm)	Part No.
DC-Red	160x110	G381321030
DC-Red	50x50	G381321036
DC-Red	25x25	G381321033
DC-Red	Ø=50	G381321035
DC-Red	Ø=25	G381321032
DC-Green	160x110	G381322030
DC-Green	50x50	G381322036
DC-Green	25x25	G381322033
DC-Green	Ø=50	G381322035
DC-Green	Ø=25	G381322032
DC-Blue	160x110	G381323030
DC-Blue	50x50	G381323036
DC-Blue	25x25	G381323033
DC-Blue	Ø=50	G381323035
DC-Blue	Ø=25	G381323032

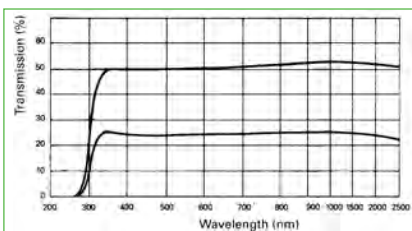Typical transmittance curve of neutral-density filters

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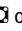
Neutral-Density Filters

Attenuate polychromatic light without altering its spectral distribution
 Transmittance $T = \Phi_T / \Phi_0$
 Optical density $D = \log 1 / T$

- Metallic films vacuum-evaporated onto glass substrates
- Hard
- Durable
- Firmly adhering
- Mounted and unmounted versions available
- Mounted types fit to Microbench  or Nanobench 
- Thickness of all filters: 2 mm

Neutralfilter
Neutral-density filters

Fassung	Ø (mm)	Transmission T (%)	T	Optische Dichte D	Toleranz Optische Dichte	Part No.
<i>Mount</i>	<i>Ø (mm)</i>	<i>Transmission T (%)</i>	<i>T</i>	<i>Optical density D</i>	<i>Tolerance optical density</i>	<i>Part No.</i>
-	12.7	50	0.5	0.3	±0.03	G371190000
-	12.7	40	0.4	0.4	±0.04	G371198000
-	12.7	20	0.2	0.7	±0.07	G371191000
-	12.7	10	0.1	1.0	±0.1	G371192000
-	12.7	5	0.05	1.3	±0.1	G371193000
-	12.7	2	0.02	1.7	±0.2	G371194000
-	12.7	1	0.01	2.0	±0.2	G371195000
-	12.7	0.1	0.001	3.0	±0.5	G371196000
-	22.4	50	0.5	0.3	±0.03	G371140000
-	22.4	40	0.4	0.4	±0.04	G371148000
-	22.4	20	0.2	0.7	±0.07	G371141000
-	22.4	10	0.1	1.0	±0.1	G371142000
-	22.4	5	0.05	1.3	±0.1	G371143000
-	22.4	2	0.02	1.7	±0.2	G371144000
-	22.4	1	0.01	2.0	±0.2	G371145000
-	22.4	0.1	0.001	3.0	±0.5	G371146000
-	50	50	0.5	0.3	±0.03	G371150000
-	50	40	0.4	0.4	±0.04	G371158000
-	50	20	0.2	0.7	±0.07	G371151000
-	50	10	0.1	1.0	±0.1	G371152000
-	50	5	0.05	1.3	±0.1	G371153000
-	50	2	0.02	1.7	±0.2	G371154000
-	50	1	0.01	2.0	±0.2	G371155000
-	50	0.1	0.001	3.0	±0.5	G371156000
☒ Mount N, clear Ø 11.5 mm	12.7	50	0.5	0.3	±0.03	G052501000
☒ Mount N, clear Ø 11.5 mm	12.7	40	0.4	0.4	±0.04	G052509000
☒ Mount N, clear Ø 11.5 mm	12.7	20	0.2	0.7	±0.07	G052502000
☒ Mount N, clear Ø 11.5 mm	12.7	10	0.1	1.0	±0.1	G052503000
☒ Mount N, clear Ø 11.5 mm	12.7	5	0.05	1.3	±0.1	G052504000
☒ Mount N, clear Ø 11.5 mm	12.7	2	0.02	1.7	±0.2	G052505000
☒ Mount N, clear Ø 11.5 mm	12.7	1	0.01	2.0	±0.2	G052506000
☒ Mount N, clear Ø 11.5 mm	12.7	0.1	0.001	3.0	±0.5	G052507000
☒ Mount CL, clear Ø 21.4 mm	22.4	50	0.5	0.3	±0.03	G063460000
☒ Mount CL, clear Ø 21.4 mm	22.4	40	0.4	0.4	±0.04	G063468000
☒ Mount CL, clear Ø 21.4 mm	22.4	20	0.2	0.7	±0.07	G063461000
☒ Mount CL, clear Ø 21.4 mm	22.4	10	0.1	1.0	±0.1	G063462000
☒ Mount CL, clear Ø 21.4 mm	22.4	5	0.05	1.3	±0.1	G063463000
☒ Mount CL, clear Ø 21.4 mm	22.4	2	0.02	1.7	±0.2	G063464000
☒ Mount CL, clear Ø 21.4 mm	22.4	1	0.01	2.0	±0.2	G063465000
☒ Mount CL, clear Ø 21.4 mm	22.4	0.1	0.001	3.0	±0.5	G063466000

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☒ / ☒ In mount fitting directly to Nanobench / Microbench, see chapter "Mounts, Posts and Holders", section "Lens Mounts" for mount dimensions.

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
Neutral Wedges

- Fabricated from neutral-density filter glass
- Cemented onto compensating plate
- Transmittance range of 1-80 %
- Faces parallel to within 10'

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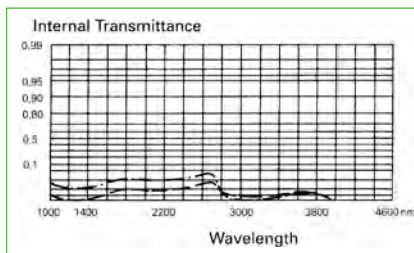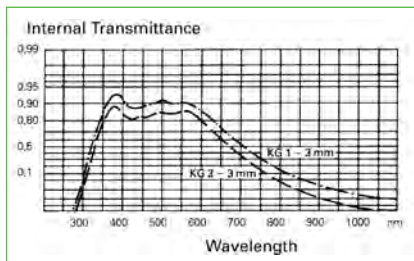
Heat-Absorbing Filters

- Fabricated from tempered glass
- Virtually colorless
- Unmounted and mounted versions available
- In mount fitting directly to Microbench

Tolerances:

- Ø 22.2mm: ± 0.2 mm; Ø 50 mm: +0/-0.6
- Thickness: ±0.2 mm
- Bubbles: 1/5x0.25
- Striations: 2/03
- Surface Finish: 5/3x0.4
- Surface Parallelism: 15'

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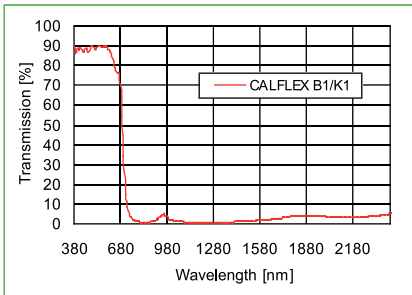
In mount fitting directly to Microbench, see chapter "Mounts, Posts and Holders", section "Lens Mounts" for mount dimensions.

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Heat-Protection Filters CALFLEX™ B1/K1

- Consists of IR-absorption glass and dielectric coating
- High suppression of IR-(Heat-)radiation
- High transmission in the visible range
- Heat-resistant up to 250 °C

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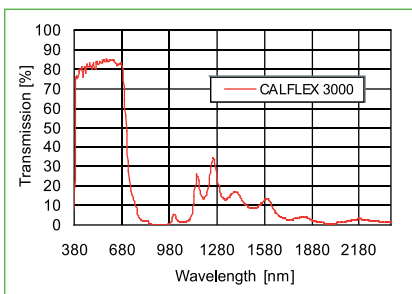
Heat-Protection Filters CALFLEX™ 3000 SP

- No absorption from dielectric coating
- Broadband heat-protection for halogen lamps
- UV-Light suppression
- High transmission in the visible range
- Heat-resistant up to 400 °C

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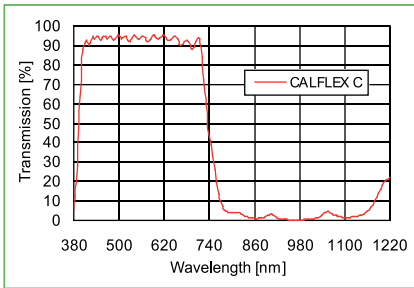
Heat-Protection Filters CALFLEX™ C


- Heat-light separation for high pressure discharge lamps
- High transmission in the visible range
- No absorption of dielectric coating
- Heat-resistant up to 400 °C

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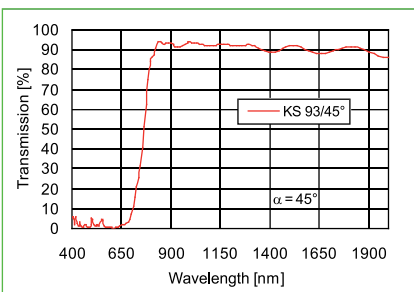Cold Light Mirror KS 93 / 45°


- For incidence angles of 45°
- Transmittance of IR light and reflection of visible light
- High reflection in the visible range
- No absorption of dielectric coating
- Heat-resistant up to 300 °C

- Thickness 1.1 mm
- $R \geq 97\%$ for 425 - 600 nm
- $T = 50\%$ for 685 ± 15 nm
- $T \geq 80\%$ for 750 - 2500 nm

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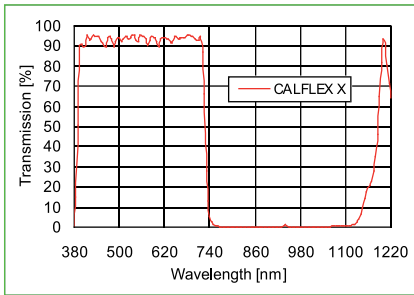
NIR-Blocking Filters CALFLEX™ X

- For photo detectors, Si-Diodes
- Very high blocking in the NIR
- High transmission in the VIS
- Heat-resistant up to 250 °C
- Thickness 1.1 mm
- $T \leq 1\%$ for 780-1100 nm
- $T = 50\%$ for 730 ±20 nm
- $T \geq 80\%$ for 420-680 nm

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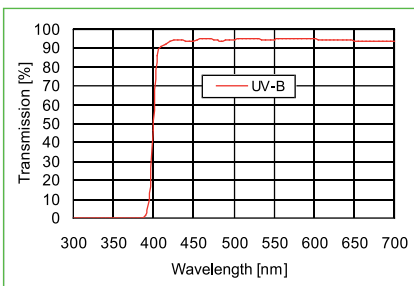UV-Blocking Filters UV-B

- Very high blocking in the UV
- High transmission in the VIS
- Heat-resistant up to 400 °C
- Thickness 1.1 mm
- $T \leq 1\%$ below 380 nm
- $T = 50\%$ for 405 ± 10 nm
- $T \geq 90\%$ average for 425 - 760 nm

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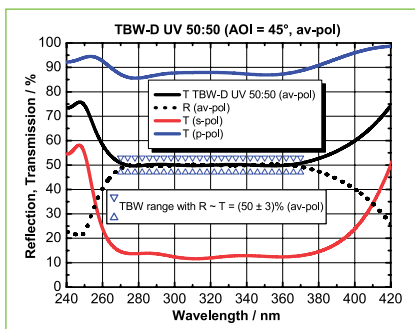
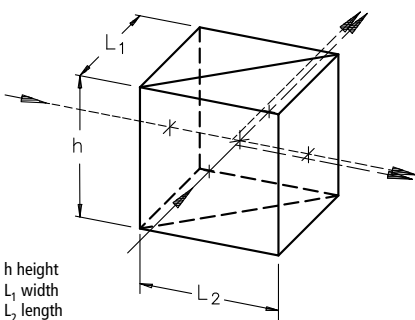
Beamsplitter Cubes UV

- For splitting or combining light beams
- No angular or lateral deflection of transmitted beams
- Reflected and transmitted beams transit identical optical paths
- Compact, rugged units
- Designed for use with low-power lasers (0.1 J / cm² for 10 ns pulses at 1064 nm)
- Fused silica
- Reflectance ≈ transmittance ±5 % for unpolarized light
- Broadband ARB1 UV coated
- 8' angular beam deflection tolerance

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

Recommended holder for cubes and prisms with 10 mm edge length:

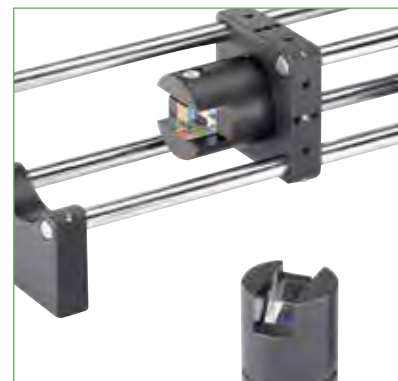Clamp holder for prism (G06 1641 000, chapter Microbench)

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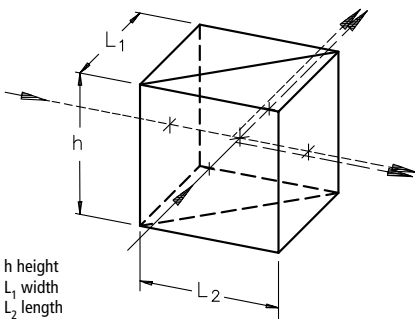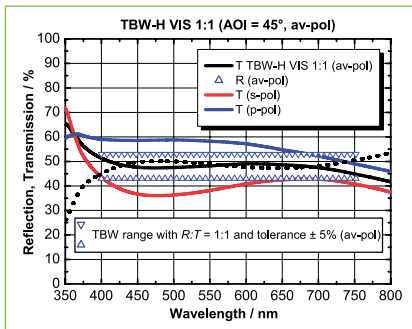
Beamsplitter Cubes VIS

- For splitting or combining light beams
- No angular or lateral deflection of transmitted beams
- Reflected and transmitted beams transit identical optical paths
- Compact, rugged units
- Designed for use with low-power lasers (0.1 J/cm² for 10 ns pulses at 1064 nm)
- N-BK7 glass
- Reflectance ≈ transmittance ± 5 % for unpolarized light
- Absorption < 5 %
- Broadband ARB2 VIS coated
- 8' angular beam deflection tolerance

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
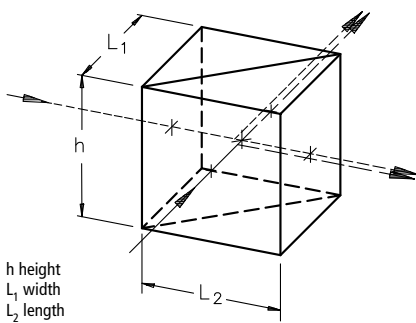
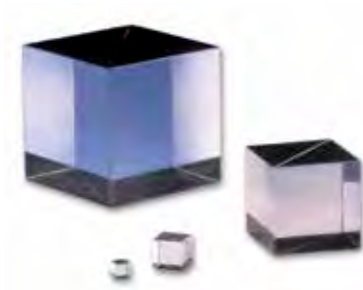
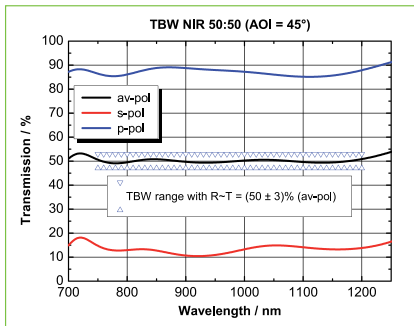
Beamsplitter Cubes NIR

- For splitting or combining light beams
- No angular or lateral deflection of transmitted beams
- Reflected and transmitted beams transit identical optical paths
- Compact, rugged units
- Designed for use with low-power lasers (0.1 J/cm² for 10 ns pulses at 1064 nm)
- N-BK7 glass
- Reflectance ≈ transmittance ±5 % for unpolarized light
- Absorption < 5 %
- Broadband AR-coated (ARB2-NIR)
- 8 arcmin angular beam deflection tolerance

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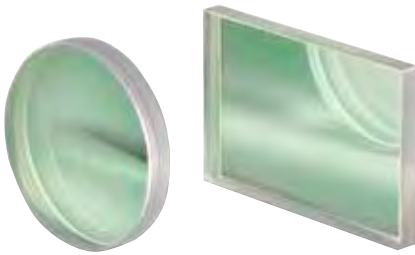
Beamsplitter Plates UV

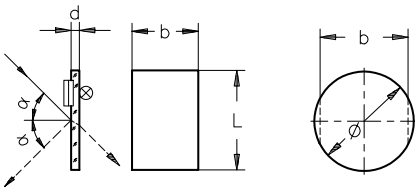
- Designed for a wavelength range of 248-400 nm
- Rectangular, circular and elliptical plates
- R : T = 20 : 80 ± 3 %
- R : T = 33 : 66 ± 4 %
- R : T = 50 : 50 ± 5 %
- AOI = 45°
- Second surface antireflective-coated with ARB2 UV

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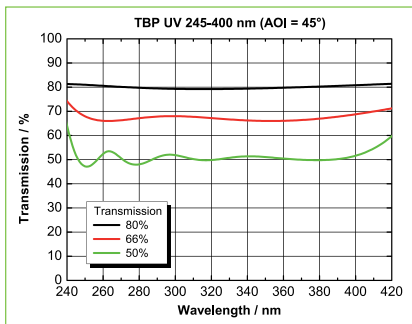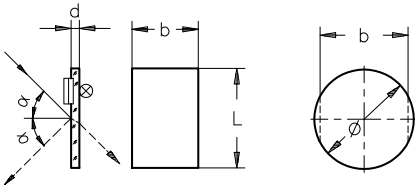
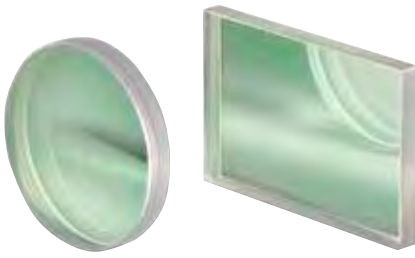
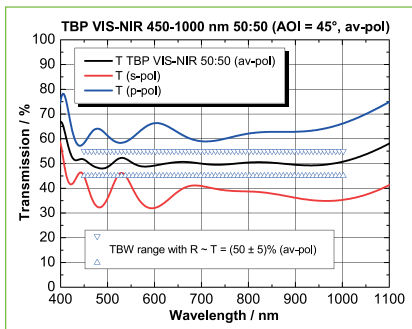
Beamsplitter Plates VIS-NIR

- For splitting or combining beams
- Also available with a wedge for eliminating etalon effects
- Second surface broadband ARB1 coated
- N-BK7 glass
- Reflectance \approx transmittance $\pm 5\%$ for unpolarized light at AOI = 45°
- Broadband, 450-1000 nm

- Other beamsplitter coatings, as well as different splitting ratios, are described in chapter Thin Film Coatings.
- For higher quantities, customized versions (other sizes, materials, coatings) available on request. Ask our sales team for a special quotation.

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Früher
In former times

Montage und Mechanikfertigung in Göttingen um 1910
Mounting and mechanics production in Goettingen about 1910

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
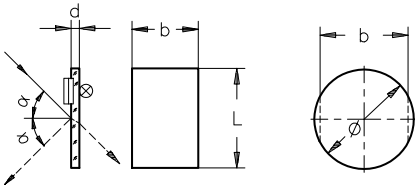
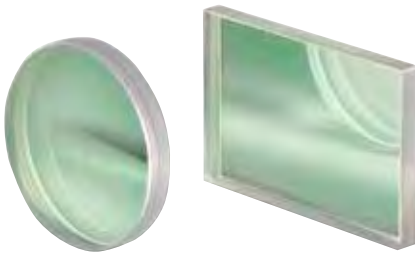
Beamsplitter Plates NIR

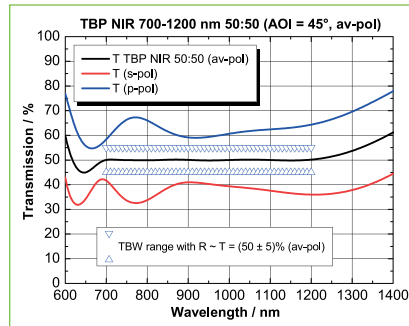- For splitting or combining beams
- Second surface broadband AR-coated (ARB2 NIR, 700-1200 nm)
- N-BK7 glass
- Reflectance \approx transmittance $\pm 5\%$ for unpolarized light at AOI = 45°
- Broadband, 700-1200 nm
- Damage Threshold: 10 J/cm² at 1064 nm, 10 ns pulse duration

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

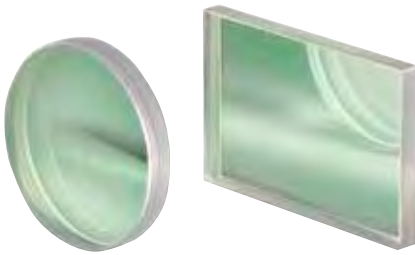
Beamsplitter Plates IR

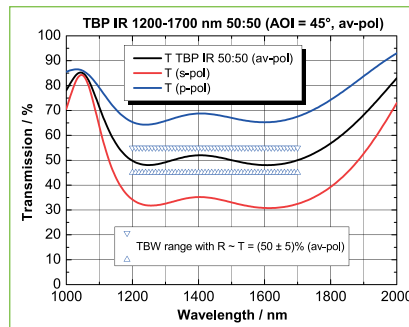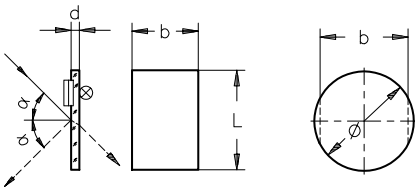
- Developed for 1250-1650 nm wavelength
- Rectangular and circular plates
- R:T = 50:50 +5 % for unpolarized light
- Second surface broadband ARB2 IR coated
- N-BK7 glass
- AOI = 45°

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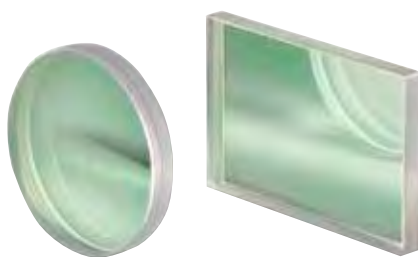
Beamsplitter Plates with Different Splitting Ratios

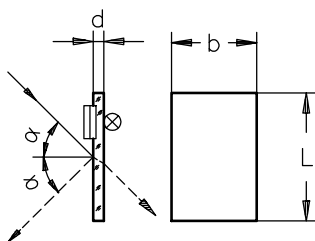- For splitting or combining beams
- Withstand high incident intensities
- Choice of different reflectance/transmittance ratios
- TBP broadband beamsplitter coating

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
- N-BK7 glass
- Second surface broadband AR-coated
- Tolerances: $b = -0.13 \text{ mm}$, $L = -0.13 \text{ mm}$, $d = \pm 0.3 \text{ mm}$
- Flatness of first surface at $633 \text{ nm } \lambda/2$
- Flatness of the back side 1λ

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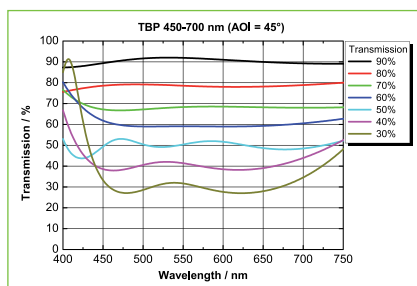
- For beam splitters 50:50, please refer to chapter "Beamsplitter Plates VIS-NIR"

Transmission ratios of Beamsplitter Plates from T=90% to T=30%, AOI = 45°

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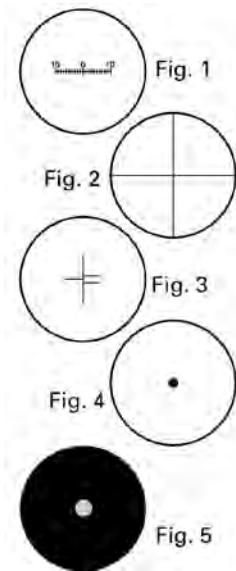
Reticles

- High-precision optical-glass rulings
- Ideally suited for positioning / measurement aids for adjusting beam paths, or for use as components in optical equipment and instruments
- Unmounted and mounted versions
- In mount fitting directly to Microbench ☒


Strichplatten (Okularmikrometer) Reticles

Fassung	Ø (mm)	Dicke d (mm)	Toleranz Ø (mm)	Teilung	Abbildung	Part No.
Mount	Ø (mm)	Thickness d (mm)	Tolerance Ø (mm)	Graduation	Fig.	Part No.
-	19	2	-0.13	5 mm in 200 divisions	1	G391122000
-	19	2	-0.13	10 mm in 200 divisions	1	G391126000
-	19	2	-0.13	dual crosshairs; 2x3 mm; 5 µm	3	G391124000
-	19	2	-0.13	crosshairs; full-width 12 µm	2	G391130000
☒	19	2	-0.13	5 mm in 200 divisions	1	G063510000
☒	19	2	-0.13	10 mm in 200 divisions	1	G063511000
☒	19	2	-0.13	crosshairs; full-width 12 µm	2	G063514000

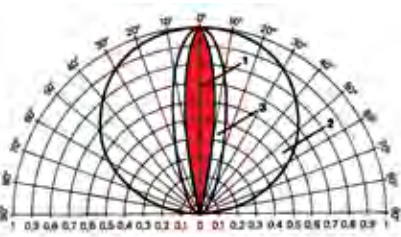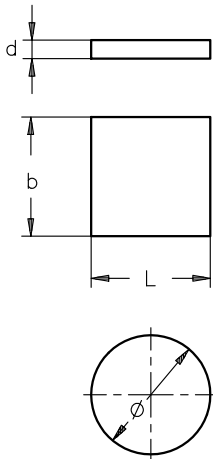Light scattering patterns:
 1. Ground glass plates,
 2. Opal glass plates

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Surface Dispersion Plates (Ground Glass)

- For observing real optical images in beam paths
- For providing uniform illumination from lighting systems
- Provide approx. 10 µm scattering center
- Other sizes available for higher quantities on request.

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Our quality criteria:

- *High-quality materials*
- *Continuous process monitoring*
- *Coatings for stringent demands in science and industry*
- *Laser-suitable coatings*
- *Expert consultation, from the development phase to the finished product*

LINOS Mirrors

Metallic, hybrid, dielectric: the standard line of Qioptiq products includes a large selection of mirrors for a variety of applications. Both broadband and highly reflective mirrors for the most common discrete wavelengths are available directly from stock. Many of the dielectric mirrors can also be used as dichroitic beam splitters.

We can prepare a quote for a tailored solution for large quantities, upon request. Just contact us!

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Special Features:

We can also supply catalog items with customized dimensions upon request.

Ideal Areas:

Laser and high-performance optics; precision optics; illumination optics; beam splitting or superimposing.

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Content

Plane Mirrors

<i>Gold Plane Mirrors, Laser Quality</i>	334
<i>Silver Rectangular Plane Mirrors, Laser Quality</i>	335
<i>Silver Rectangular Plane Mirrors, Standard Quality</i>	336
<i>Silver Circular Plane Mirrors, Laser Quality</i>	337
<i>Silver Elliptical Plane Mirrors, Laser Quality</i>	338
<i>Silver Elliptical Plane Mirrors, Standard Quality</i>	339
<i>Aluminized Plane Mirrors - Rectangular plane</i>	341
<i>Aluminized Plane Mirrors - Circular plane</i>	342
<i>Aluminized Plane Mirrors - Elliptical plane</i>	344
<i>Dielectric-Coated Plane Mirrors DLHS VUV 157 nm</i>	345
<i>Dielectric-Coated Plane Mirrors DLHS UV 248 nm</i>	346
<i>Dielectric-Coated Plane Mirrors DLHS UV 266 nm</i>	346
<i>Dielectric-Coated Plane Mirrors DLHS UV 308 nm</i>	347
<i>Dielectric-Coated Plane Mirrors DLHS UV 351-355 nm</i>	348
<i>Dielectric-Coated Plane Mirrors DLHS UV 405 nm</i>	349
<i>Dielectric-Coated Plane Mirrors DLHS 532 nm</i>	349
<i>Dielectric-Coated Plane Mirrors DLHS 633 nm</i>	350
<i>Dielectric Coated Plane Mirror DLHS 755-840 nm, Ti:Sa</i>	351
<i>Dielectric-Coated Plane Mirrors DLHS NIR 1064 nm</i>	352
<i>Dielectric-Coated Plane Mirrors DLHS NIR 1540 nm</i>	352
<i>Dielectric-Coated Plane Mirrors DLB UV 230-400 nm</i>	353
<i>Dielectric-Coated Plane Mirrors DLB SBB 350-950 nm</i>	354
<i>Dielectric-Coated Plane Mirrors LBSM-UV</i>	355
<i>Dielectric-Coated Plane Mirrors LBSM-VIS</i>	355
<i>Dielectric-Coated Plane Mirrors LBSM-NIR</i>	356
<i>Dielectric-Coated Plane Mirrors LBSM-YAG</i>	356
<i>Plane Mirror SILFLEX</i>	357

Concave Mirrors

<i>Front Surface Mirrors for Illumination Systems</i>	358
---	-----

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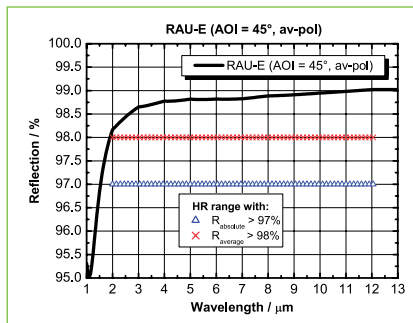
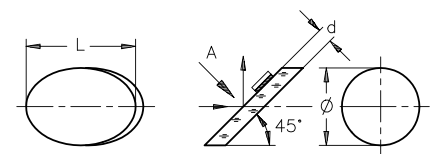
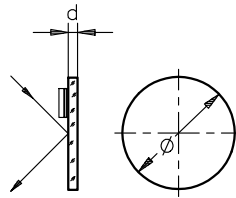
Gold Plane Mirrors, Laser Quality

- Front surface gold mirror for the IR spectral range up to 25 μm
- Resistant against environmental influences
- Constant high reflection within a broad spectral and angle-of-incidence range without absorption bands
- Average reflection > 98 % in the range of $0^\circ \leq \text{AOI} \leq 60^\circ$, unpolarized
- Low expansion material
- Dielectric protection layer for improved environmental stability

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Of course highest quality mirrors need excellent mirror mounts for high precision X-Y adjustments. Our premium class Lees mirror mounts with least possible backlash due to counter surfaces of sapphire and the special Krytox lubricant will elate you too.

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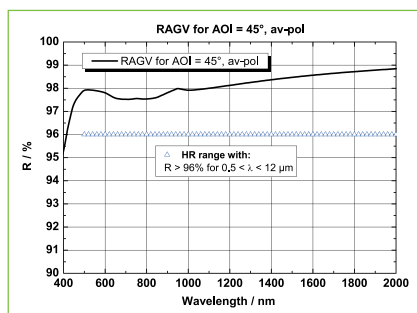
☑ Recommended mounts: Lees™ Mirror Mounts in chapter Mirror Mounts.

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
Example: RAGV (AOI=45°, unpolarized)

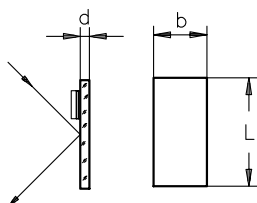
Silver Rectangular Plane Mirrors, Laser Quality

- Front surface rectangular mirrors
- Equal reflectance without absorption band
- Angle of incidence from 0° to 60°
- Resistant against environmental influences
- Double dielectric protection layer
- Suitable for short-pulse applications
- Highest reflectance for 500 nm to 12 µm wavelength range
- Guaranteed reflection ($0^\circ \leq \text{AOI} \leq 60^\circ$, unpolarized) > 96 % in the range of $500 \text{ nm} \leq \lambda \leq 780 \text{ nm}$, > 97 % in the range of $780 \text{ nm} \leq \lambda \leq 10 \mu\text{m}$
- Silver front surface coating RAGV for 500 nm up to 12 µm
- Unmounted
- Low expansion material
- Damage threshold $H_\infty > 1.5 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz

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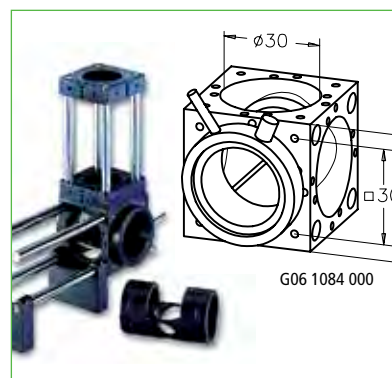Recommended mirror holder for rectangular mirrors 20x30mm for the Microbench: Cube 30 with build-in rotatable plate holder for plano optics 20 x 30 mm (Beam-deflector cube G06 1084 000, chapter Microbench).


Planspiegel, eckige Form, Laserqualität, silberbeschichtet Silver Rectangular Plane Mirrors, Laser Quality


Größe bxL (mm)	Dicke d (mm)	Toleranz d (mm)	Toleranz b (mm)	Toleranz L (mm)	Planität (λ)	Part No.
Size bxL (mm)	Thickness d (mm)	Tolerance d (mm)	Tolerance b (mm)	Tolerance L (mm)	Flatness (λ)	Part No.
10x15	2	±0.2	-0.09	-0.11	1/10	G340331400
20x30	4	±0.2	-0.13	-0.13	1/10	G340536000
60x85	8	±0.2	-0.19	-0.22	1/10	G340332400

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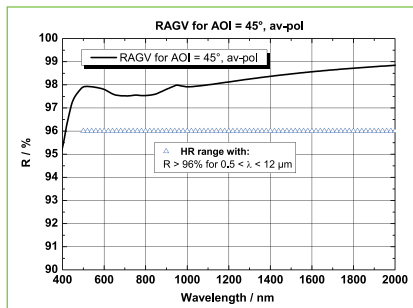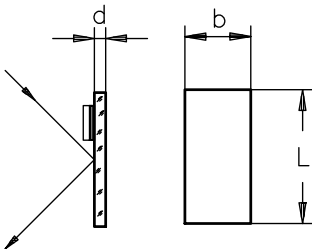Silver Rectangular Plane Mirrors, Standard Quality

- Front surface rectangular mirrors
 - Equal reflectance without absorption band
 - Angle of incidence from 0° to 60°
 - Resistant against environmental influences
 - Double dielectric protection layer
 - Suitable for short-pulse applications
- Highest reflectance for 500 nm to 12 μm wavelength range
 - Guaranteed reflection (0° ≤ AOI ≤ 60°, unpolarized) > 96 % in the range of 500 nm ≤ λ ≤ 780 nm, > 97 % in the range of 780 nm ≤ λ ≤ 10 μm
 - Damage threshold $H_{10} > 1.5 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
 - Silver front surface coating RAGV for 500 nm up to 12 μm
 - Unmounted
 - Low expansion material

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Please find further information about application examples, guidelines, frequently asked questions and more in our Brochure "Microbench, Bench and Rail systems". You will find the brochure for download at each Microbench product under "Docs + Drawings".

Planspiegel, eckige Form, Standardqualität, silberbeschichtet Silver Rectangular Plane Mirrors, Standard Quality

Größe bxL (mm)	Dicke d (mm)	Toleranz d (mm)	Toleranz b (mm)	Toleranz L (mm)	Planität (λ)	Part No.
Size bxL (mm)	Thickness d (mm)	Tolerance d (mm)	Tolerance b (mm)	Tolerance L (mm)	Flatness (λ)	Part No.
10x10	2	±0.2	-0.09	-0.09	1/4	G340611000
10x15	2	±0.2	-0.09	-0.11	1/2	G340614000
15x20	2	±0.2	-0.11	-0.13	1/2	G340615000
20x30	2.5	±0.2	-0.13	-0.13	1/2	G340636000
40x58	8	±0.2	-0.16	-0.19	1/4	G340642000
60x85	8	±0.2	-0.19	-0.22	1	G340662000
80x120	8	±0.2	-0.19	-0.22	2	G340682000
100x150	10	±0.2	-0.22	-0.25	4	G34066400

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
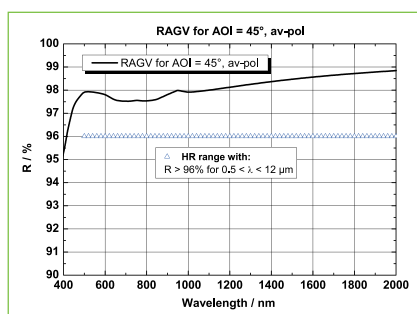
Silver Circular Plane Mirrors, Laser Quality

- Front surface rectangular mirrors
- Equal reflectance without absorption band
- Angle of incidence from 0° to 60°
- Resistant against environmental influences
- Double dielectric protection layer
- Suitable for short-pulse applications

- Highest reflectance for 500 nm to 12 µm wavelength range
- Guaranteed reflection ($0^\circ \leq \text{AOI} \leq 60^\circ$, unpolarized) > 96 % in the range of $500 \text{ nm} \leq \lambda \leq 780 \text{ nm}$, > 97 % in the range of $780 \text{ nm} \leq \lambda \leq 10 \mu\text{m}$
- Damage threshold $H_{10} > 1.5 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Silver front surface coating RAGV for 500 nm up to 12 µm
- Unmounted
- Low expansion material

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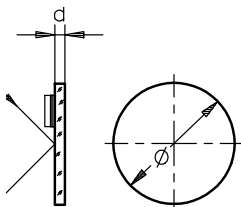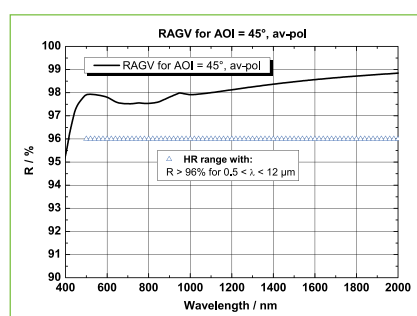
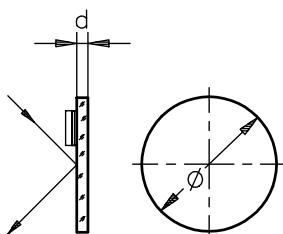
Silver Circular Plane Mirrors, Standard Quality

- Front surface rectangular mirrors
- Equal reflectance without absorption band
- Angle of incidence from 0° to 60°
- Resistant against environmental influences
- Double dielectric protection layer
- Suitable for short-pulse applications

- Highest reflectance for 500 nm to 12 μm wavelength range
- Guaranteed reflection ($0^\circ \leq \text{AOI} \leq 60^\circ$, unpolarized) > 96 % in the range of $500 \text{ nm} \leq \lambda \leq 780 \text{ nm}$, > 97 % in the range of $780 \text{ nm} \leq \lambda \leq 10 \mu\text{m}$
- Damage threshold $H_{90} > 1.5 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Silver front surface coating RAGV for 500 nm up to 12 μm
- Unmounted
- Low expansion material

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Qioptiq uses glass from established manufacturers in the production of precision optics components. This ensures the consistent high quality of our lenses, plano-optics and mirrors.

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☒ Recommended mounts: Lees™ Mirror Mounts in chapter Mirror Mounts.

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
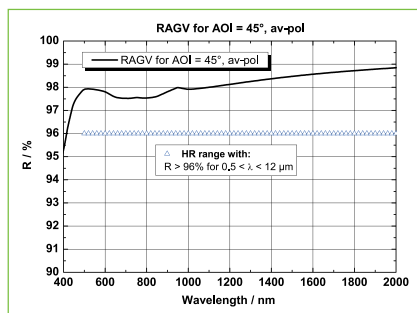
Silver Elliptical Plane Mirrors, Laser Quality

- Front surface rectangular mirrors
- Equal reflectance without absorption band
- Angle of incidence from 0° to 60°
- Resistant against environmental influences
- Double dielectric protection layer
- Suitable for short-pulse applications
- Highest reflectance for 500 nm to 12 µm wavelength range
- Guaranteed reflection ($0^\circ \leq \text{AOI} \leq 60^\circ$, unpolarized) > 96 % in the range of $500 \text{ nm} \leq \lambda \leq 780 \text{ nm}$, > 97 % in the range of $780 \text{ nm} \leq \lambda \leq 10 \mu\text{m}$
- Silver front surface coating RAGV for 500 nm up to 12 µm
- Unmounted
- Low expansion material
- Damage threshold $H_{\infty} > 1.5 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz

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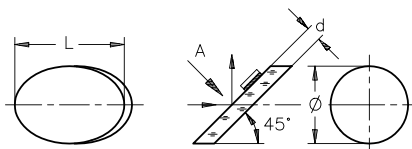
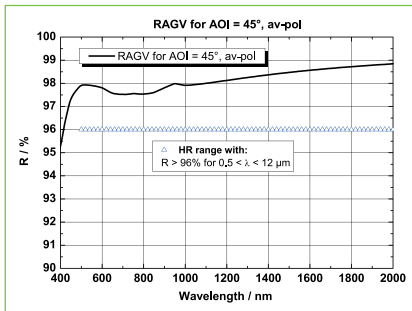
Example: RAGV (AOI=45°, unpolarized)

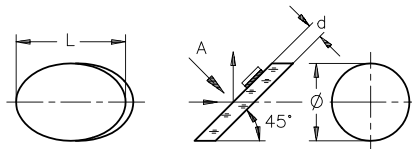Silver Elliptical Plane Mirrors, Standard Quality

- Front surface rectangular mirrors
- Equal reflectance without absorption band
- Angle of incidence from 0° to 60°
- Resistant against environmental influences
- Double dielectric protection layer Suitable for short-pulse applications

- Highest reflectance for 500 nm to 12 μm wavelength range
- Guaranteed reflection ($0^\circ \leq \text{AOI} \leq 60^\circ$, unpolarized) $> 96\%$ in the range of $500 \text{ nm} \leq \lambda \leq 780 \text{ nm}$, $> 97\%$ in the range of $780 \text{ nm} \leq \lambda \leq 10 \mu\text{m}$
- Damage threshold $H_{\infty} > 1.5 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Silver front surface coating RAGV for 500 nm up to 12 μm
- Unmounted
- Low expansion material


Planspiegel, ovale Form, Standardqualität, silberbeschichtet Silver Elliptical Plane Mirrors, Standard Quality

Größe ØxL (mm)	Dicke d (mm)	Toleranz d (mm)	Toleranz Ø (mm)	Planität (λ)	Part No.
Size ØxL (mm)	Thickness d (mm)	Tolerance d (mm)	Tolerance Ø (mm)	Flatness (λ)	Part No.
10x14	2	± 0.2	-0.06	1/2	G340612000
12.5x18	2	± 0.2	-0.08	1/2	G340616000
22.4x31.5	3.5	± 0.2	-0.10	1/2	G340623000
31.5x45	4	± 0.2	-0.12	1/2	G340632000
50x71	6	± 0.2	-0.12	1	G340651000
80x113	12.5	± 0.2	-0.14	1/4	G340681000

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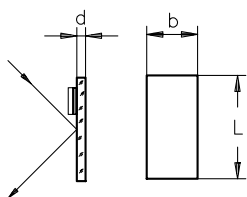The substrate low expansion material is a glass with $n_d = 1.47$ and a coefficient of expansion $\alpha_{20-300 \text{ }^\circ\text{C}} = 3.3 \cdot 10^{-6} / \text{K}$ (i.e. Pyrex, DURAN, Tempax, Borofloat).

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

Aluminized Plane Mirrors - Rectangular plane

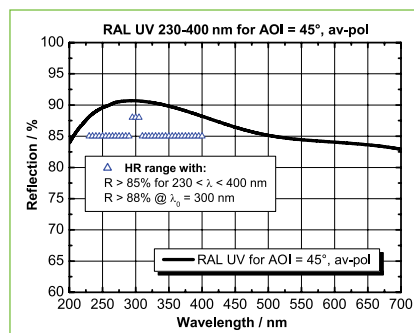- Standard aluminium mirrors for the spectral ranges VIS to IR up to $10 \mu\text{m}$ or UV to $3 \mu\text{m}$, respectively
- High reflection within a broad spectral and angle-of-incidence range
- Protected by an SiO_2 overcoating (RAL) or MgF_2 protective overcoating for the UV-region (RAL UV)
- Low expansion material
- RAL-UV: Guaranteed reflection $> 85\%$ in the range of $230 \text{ nm} \leq \lambda \leq 400 \text{ nm}$, $> 88\%$ at $\lambda_0 = 300 \text{ nm}$ for $\text{AOI} = 45^\circ$, unpolarized
Damage threshold $H_{\infty} > 200 \text{ mJ/cm}^2$ at 193 nm with 30 ns laser pulses (s-on-1) at 10 Hz
- RAL VIS: Guaranteed reflection $> 85\%$ in the range of $400 \text{ nm} \leq \lambda \leq 700 \text{ nm}$, $> 88\%$ at $\lambda_0 = 550 \text{ nm}$ for $\text{AOI} = 45^\circ$, unpolarized
Damage threshold $H_{\infty} > 1.5 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz

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
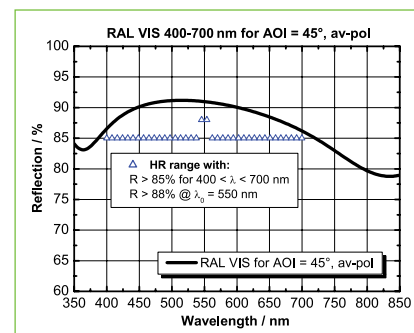
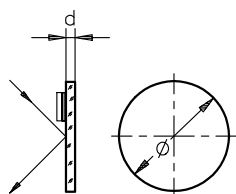
Aluminized Plane Mirrors - Circular plane

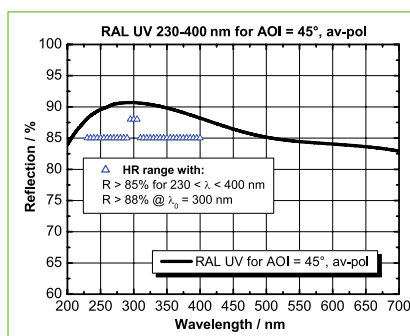- Standard aluminium mirrors for the spectral ranges VIS to IR up to $10 \mu\text{m}$ or UV to $3 \mu\text{m}$, respectively
- High reflection within a broad spectral and angle-of-incidence range
- Protected by an SiO_2 overcoating (RAL) or MgF_2 protective overcoating for the UV-region (RAL UV)
- Low expansion material
- RAL-UV: Guaranteed reflection $> 85\%$ in the range of $230 \text{ nm} \leq \lambda \leq 400 \text{ nm}$, $> 88\%$ at $\lambda_0 = 300 \text{ nm}$ for $\text{AOI} = 45^\circ$, unpolarized
Damage threshold $H_{\infty} > 200 \text{ mJ/cm}^2$ at 193 nm with 30 ns laser pulses (s-on-1) at 10 Hz
- RAL VIS: Guaranteed reflection $> 85\%$ in the range of $400 \text{ nm} \leq \lambda \leq 700 \text{ nm}$, $> 88\%$ at $\lambda_0 = 550 \text{ nm}$ for $\text{AOI} = 45^\circ$, unpolarized
Damage threshold $H_{\infty} > 1.5 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz

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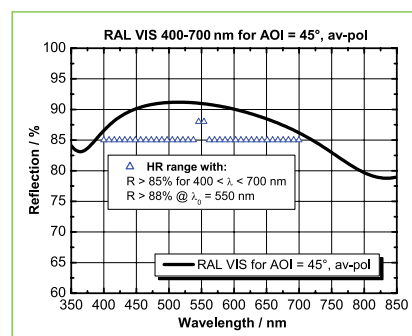The mounts and adjustment units you need for mounting mirrors in our bench systems are listed in chapter "Mounts and Posts", "Microbench", and "Mirror Mounts":

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

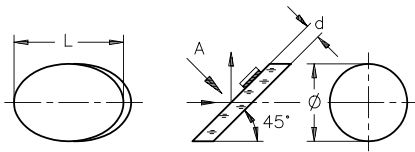
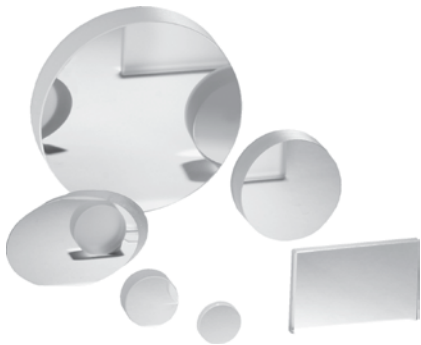
Aluminized Plane Mirrors - Elliptical plane

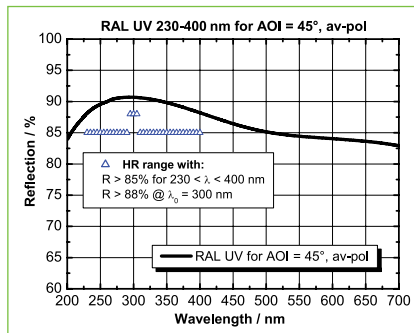- Standard aluminium mirrors for the spectral ranges VIS to IR up to $10 \mu\text{m}$ or UV to $3 \mu\text{m}$, respectively
- High reflection within a broad spectral and angle-of-incidence range
- Protected by an SiO_2 overcoating (RAL) or MgF_2 protective overcoating for the UV-region (RAL UV)
- Low expansion material
- RAL-UV: Guaranteed reflection $> 85\%$ in the range of $230 \text{ nm} \leq \lambda \leq 400 \text{ nm}$, $> 88\%$ at $\lambda_0 = 300 \text{ nm}$ for $\text{AOI} = 45^\circ$, unpolarized
Damage threshold $H_{\infty} > 200 \text{ mJ/cm}^2$ at 193 nm with 30 ns laser pulses (s-on-1) at 10 Hz
- RAL VIS: Guaranteed reflection $> 85\%$ in the range of $400 \text{ nm} \leq \lambda \leq 700 \text{ nm}$, $> 88\%$ at $\lambda_0 = 550 \text{ nm}$ for $\text{AOI} = 45^\circ$, unpolarized
Damage threshold $H_{\infty} > 1.5 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz

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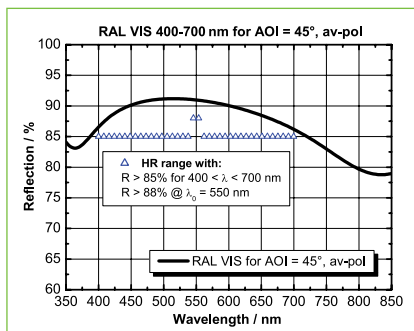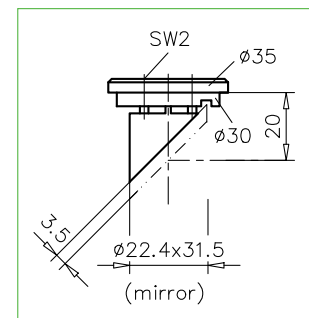
G06 3730 000

Recommended mirror holders for elliptical mirrors 22.4 x 31.5 mm:

- Adjustable mirror insert 30 (G 06 3730 000) to mount in Cube 30 of Microbench
- Beamsteering mirror mount, adjustable (G06 5096 000) to integrate in Beam deflector cube C40 of our Tube Mounting System or in Microbench mounting plates 30

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
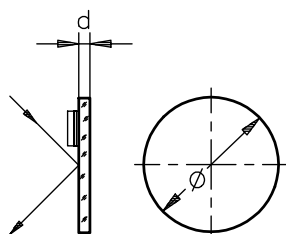
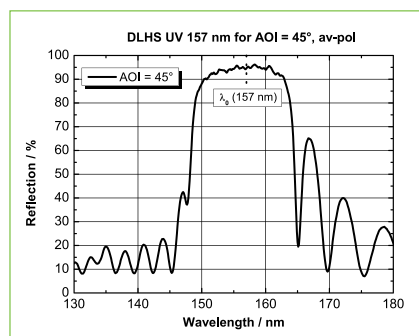
Dielectric-Coated Plane Mirrors DLHS VUV 157 nm

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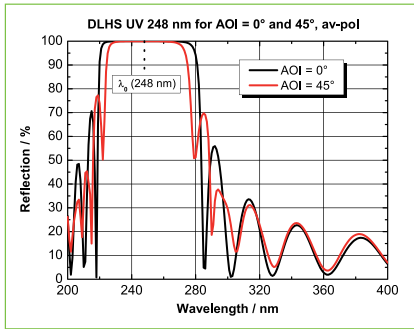DLHS DUV for 248 nm (AOI = 0°) and DLHS DUV for 248 nm (AOI = 45°), unpolarized

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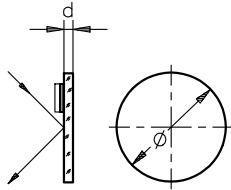
Dielectric-Coated Plane Mirrors DLHS UV 248 nm


- For high-power / high-energy laser applications
- Mirror with dielectric UV High-power coating
- Beam steering mirror for 45° angle of incidence, 90° beam deflection
- Polished back surface
- Extreme low scattering losses

- Flatness $\lambda/10$
- Fused silica substrate
- Front surface RMS ≤ 1 nm polished
- Guaranteed reflection $> 99\%$ at 248 nm, both with AOI = 0° or AOI = 45° (average polarization)
- $> 80\%$ transmission for $\lambda > 300$ nm (additional AR coating required)
- Damage thresholds LIDT > 1 J/cm² at 248 nm with 30 ns laser pulses 1000-on-1

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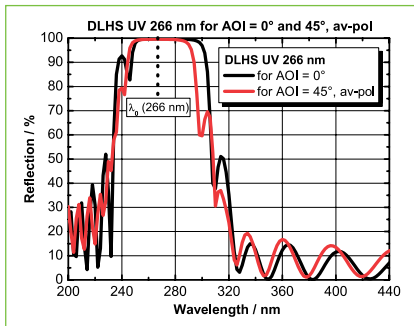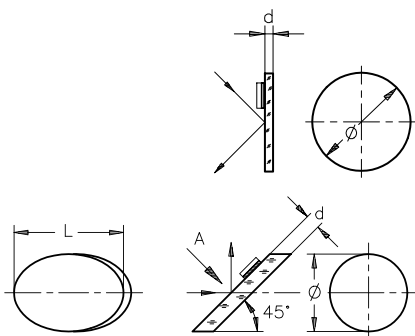
Dielectric-Coated Plane Mirrors DLHS UV 266 nm


- For high-power / high-energy laser applications
- Mirror with dielectric VUV High-power coating
- Also suitable as dichroic / longpass
- Beam steering mirror for 45° angle of incidence, 90° beam deflection
- Polished back surface
- Extreme low scattering losses

- Flatness $\lambda/10$
- Fused silica substrate
- Front surface RMS ≤ 1 nm polished
- Guaranteed reflection $> 99.5\%$ at 266 nm, both with AOI = 0° or AOI = 45° (average polarization)
- $> 80\%$ transmission for $\lambda > 300$ nm (additional AR coating required)
- Damage thresholds $H_{\infty} > 5$ J/cm² at 308 nm with 15 ns laser pulses (s-on-1) at 10 Hz

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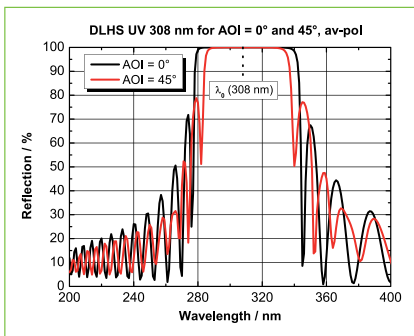
DLHS UV for 308 nm (AOI = 0°) and DLHS UV for 308 nm (AOI = 45°), unpolarized

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Dielectric-Coated Plane Mirrors DLHS UV 308 nm

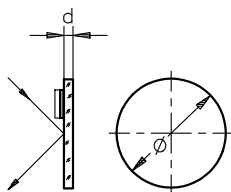- For high-power / high-energy laser applications
- Mirror with dielectric VUV High-power coating
- Also suitable as dichroic / longpass
- Beam steering mirror for 45° angle of incidence, 90° beam deflection
- Polished back surface
- Extreme low scattering losses
- Flatness $\lambda/10$
- Fused silica substrate
- Front surface $RMS \leq 1$ nm polished
- Guaranteed reflection $> 99.5\%$ at 308 nm, both with AOI = 0° or AOI = 45° (average polarization)
- $>80\%$ transmission for $\lambda > 400$ nm (additional AR coating required)
- Damage thresholds $H_{10} > 5$ J/cm² at 308 nm with 15 ns laser pulses (s-on-1) at 10 Hz

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
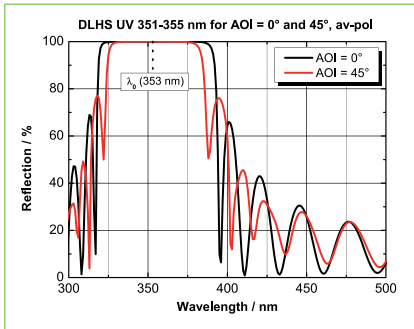
Please find further information about application examples, guidelines, frequently asked questions and more in our Brochure "Microbench, Bench and Rail systems. You will find the brochure for download at each Microbench product under "Docs + Drawings".

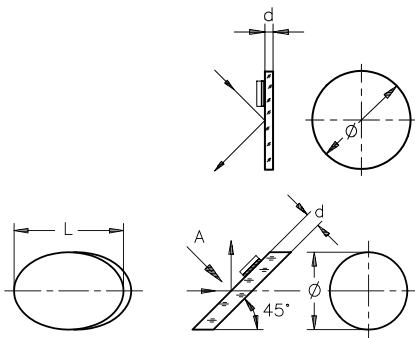DLHS UV for 351-355 nm (AOI = 0°) and DLHS UV for 351-355 nm (AOI = 45°), unpolarized

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Dielectric-Coated Plane Mirrors DLHS UV 351-355 nm

- Mirror with dielectric VUV High-power coating
- For high-power/high-energy laser applications
- Beam steering mirror for 45° angle of incidence, 90° beam deflection
- Also suitable as dichroic / longpass
- Polished back surface
- Extreme low scattering losses

- Flatness $\lambda/10$
- Fused silica substrate
- Front surface $RMS \leq 1 \text{ nm polished}$
- Guaranteed reflection $> 99.5 \%$ at 351-355 nm, both with AOI = 0° or AOI = 45° (average polarization)
- $> 80 \%$ transmission for $\lambda > 450 \text{ nm}$ (additional AR coating required)
- Damage thresholds $H_{\infty} > 5 \text{ J/cm}^2$ at 308 nm with 15 ns laser pulses (s-on-1) at 10 Hz

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*) Recommended mirror mounts for elliptical mirrors 22.4x31.5 mm:
 - Adjustable mirror insert 30 (G 06 3730 000) to mount in Cube 30, described in chapter Microbench
 - Beamsteering mirror mount, adjustable (G06 5096 000) described in chapter Tube Mounting System

Selection guide

Passende Fassungen und Justierungen zur Montage von Spiegeln finden Sie in den Kapiteln Aufbaumaterial, Mikrobank und Spiegelhalter:
The mounts and adjustment units you need for mounting mirrors in our bench systems are listed in chapter "Mounts and Posts", "Microbench", and "Mirror Mounts":

G03 6185 000

G03 6058 000

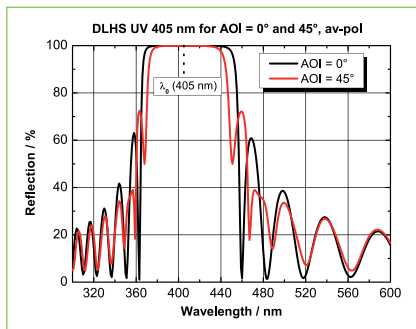
G06 5002 000

G03 5916 000


Lees Spiegelhalter
Lees Mirror Mounts

adjust.x Spiegelhalter
adjust.x Mirror Mounts

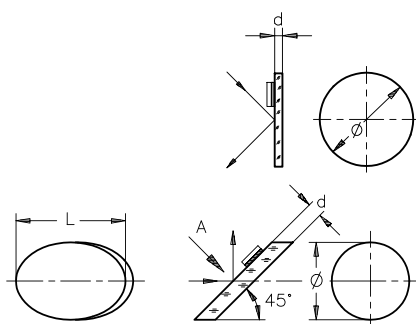DLHS UV for 405 nm (AOI = 0°) and DLHS UV for 405 nm (AOI = 45°), unpolarized

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
Dielectric-Coated Plane Mirrors DLHS UV 405 nm

- Mirror with dielectric UV High-power coating
- For high-power/high-energy laser applications
- Beam steering mirror for 0 - 50° angle of incidence
- Also suitable as dichroit / longpass
- Polished back surface
- Extreme low scattering losses


- Flatness $\lambda/10$
- Fused silica substrate
- Front surface $RMS \leq 1 \text{ nm}$ polished
- Guaranteed reflection $> 99.5 \%$ at 405 nm, both with AOI = 0° or AOI = 45° (average polarization)
- Damage thresholds $H_{\infty} > 5 \text{ J/cm}^2$ at 308 nm with 15 ns laser pulses (s-on-1) at 10 Hz

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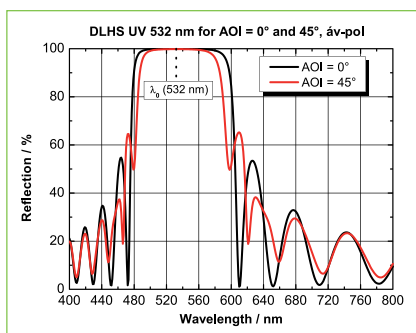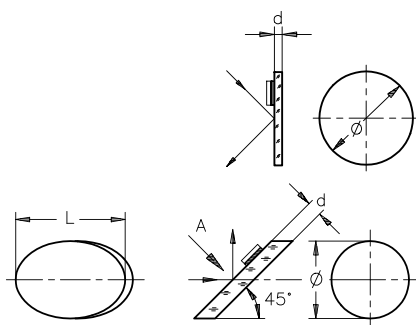
Dielectric-Coated Plane Mirrors DLHS 532 nm

- For high-power / high-energy laser applications
- Mirror with dielectric high power coating
- Beam steering mirror for 45° angle of incidence, 90° beam deflection
- Also suitable as dichroit / longpass
- Polished back surface
- Extreme low scattering losses


- Flatness $\lambda/10$
- Fused silica substrate
- Front surface $RMS \leq 1 \text{ nm}$ polished
- Guaranteed reflection $> 99.8 \%$ for 532 nm for AOI = 0° or AOI = 45° (average polarization)
- $> 90 \%$ transmission for $\lambda > 750 \text{ nm}$ (additional AR coating required)
- Damage threshold $H_{\infty} > 10 \text{ J/cm}^2$ at 532 nm with 15 ns laser pulses (s-on-1) at 10 Hz

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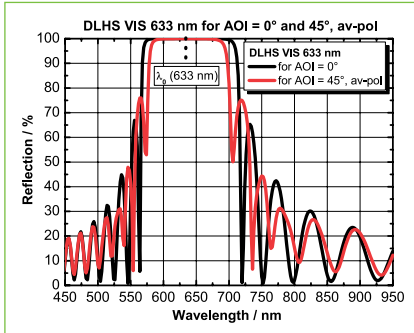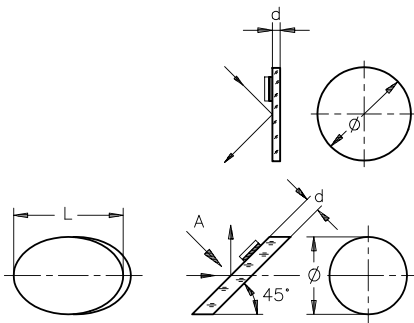Dielectric-Coated Plane Mirrors DLHS 633 nm

- For high-power / high-energy laser applications
- Mirror with dielectric high power coating
- Beam steering mirror for 45° angle of incidence, 90° beam deflection
- Also suitable as Dichroic / longpass
- Polished back surface
- Extreme low scattering losses

- Flatness $\lambda/10$
- Fused silica substrate
- Front surface $RMS \leq 1$ nm polished
- Guaranteed reflection $> 99.8\%$ for 633 nm for AOI = 0° or AOI = 45° (average polarization)
- $>90\%$ transmission for $\lambda > 850$ nm (additional AR coating required)
- Damage threshold $H_{\infty} > 10$ J / cm² at 532 nm with 15 ns laser pulses (s-on-1) at 10 Hz

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*) Recommended mirror mounts for elliptical mirrors 22.4x31.5 mm:
- Adjustable mirror insert 30 (G 06 3730 000) to mount in Cube 30, described in chapter Microbench
- Beamsteering mirror mount, adjustable (G06 5096 000) described in chapter Tube Mounting System

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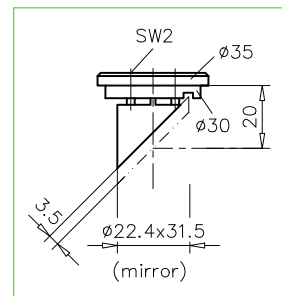
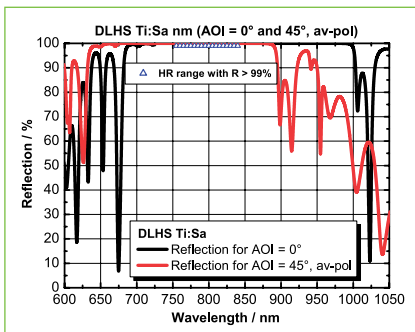
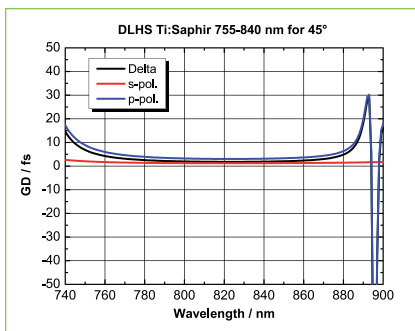
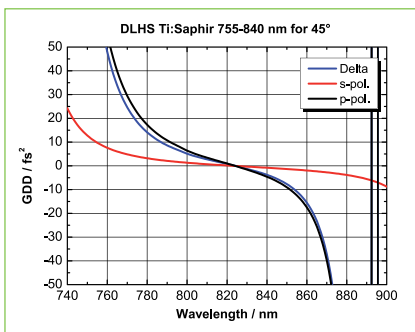
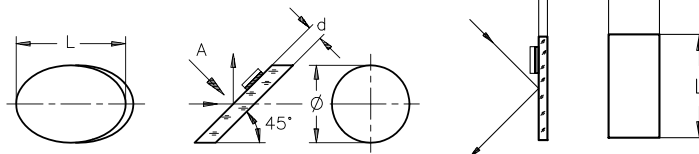
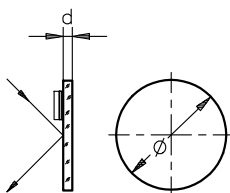
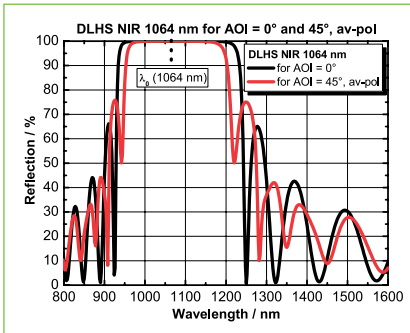
*) Recommended mirror mounts for elliptical mirrors 22.4x31.5 mm:
 - Adjustable mirror insert 30 (G 06 3730 000) to mount in Cube 30, described in chapter Microbench
 - Beamsteering mirror mount, adjustable (G06 5096 000) described in chapter Tube Mounting System


Dielectric Coated Plane Mirror DLHS 755-840 nm, Ti:Sa

- For high-power/high-energy applications
- With broadband dielectric coating designed for Ti:Sapphire lasers
- Polished back surface
- Extreme low scattering losses
- Use at 0-45° angle of incidence

- Flatness $\lambda/10$
- Fused silica substrates
- Guaranteed reflection > 99 % for the HR range 755 nm to 840 nm for $0^\circ \leq AOI \leq 45^\circ$ (average polarization)
- Damage threshold $H_{10} > 10 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz

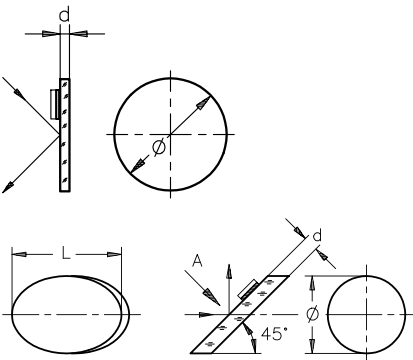DLHS NIR for 1064 nm (AOI = 0°) and DLHS NIR for 1064 nm (AOI = 45°), unpolarized

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Dielectric-Coated Plane Mirrors DLHS NIR 1064 nm

- For high-power / high-energy laser applications
- Mirror with dielectric high-power NIR coating
- Beam steering mirror for 45° angle of incidence, 90° beam deflection
- Also suitable as dichroic / shortpass
- Polished back surface
- Extreme low scattering losses

- Flatness $\lambda/10$
- Fused silica substrate
- Front surface $RMS \leq 1\text{ nm}$ polished
- Guaranteed reflection $> 99.8\%$ for 1064 nm for AOI = 0° or AOI = 45° (average polarization)
- Damage threshold $H_{\infty} > 30\text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz

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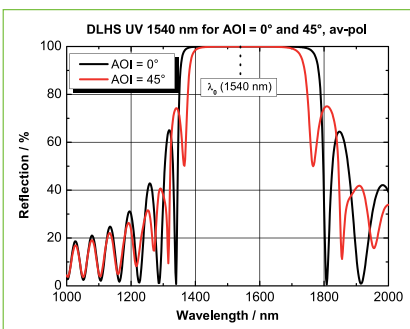
*) Recommended mirror mounts for elliptical mirrors 22.4x31.5 mm:
 - Adjustable mirror insert 30 (G 06 3730 000) to mount in Cube 30, described in chapter Microbench
 - Beamsteering mirror mount, adjustable (G06 5096 000) described in chapter Tube Mounting System

DLHS NIR for 1540 nm (AOI = 0°) and DLHS NIR for 1540 nm (AOI = 45°), unpolarized

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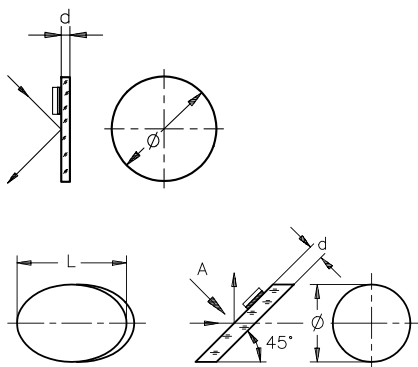
Dielectric-Coated Plane Mirrors DLHS NIR 1540 nm

- For high-power / high-energy laser applications
- Mirror with dielectric High-power NIR coating
- Beam steering mirror for 45° angle of incidence, 90° beam deflection
- Also suitable as dichroic / shortpass
- Polished back surface
- Extreme low scattering losses


- Flatness $\lambda/10$
- Fused silica substrate
- Front surface $RMS \leq 1\text{ nm}$ polished
- Guaranteed reflection $> 99.8\%$ for 1540 nm for AOI = 0° or AOI = 45° (average polarization)
- Damage threshold $H_{\infty} > 30\text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz

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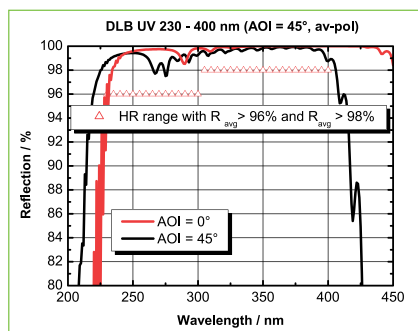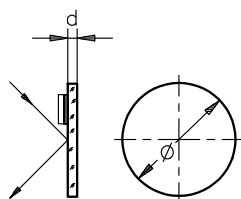
Dielectric-Coated Plane Mirrors DLB UV 230-400 nm

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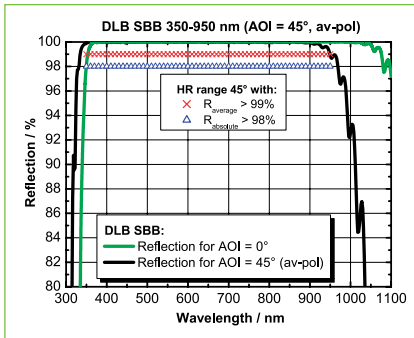
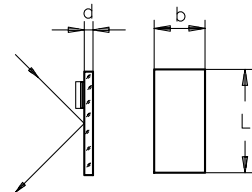
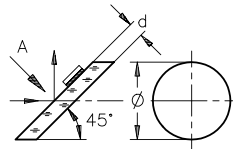
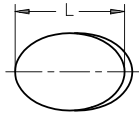
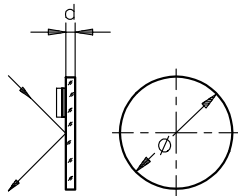
Dielectric-Coated Plane Mirrors DLB SBB 350-950 nm

- For high-power / high-energy laser applications
- Dielectric laser mirror
- Polished back surface (only fused silica)
- Extreme low scattering losses
- Beam steering mirror for 45° angle of incidence, 90° beam deflection

- Fused silica substrate
- R abs > 98 % at 350-950 nm (AOI = 45°, unpolarized) and at 400-1050 nm (AOI = 0°, unpolarized)
- R avg > 99 % at 350-950 nm (AOI = 45°, unpolarized) and at 400-1050 nm (AOI = 0°, unpolarized)

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

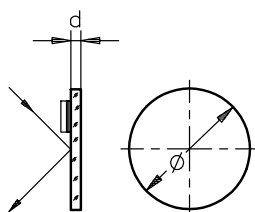
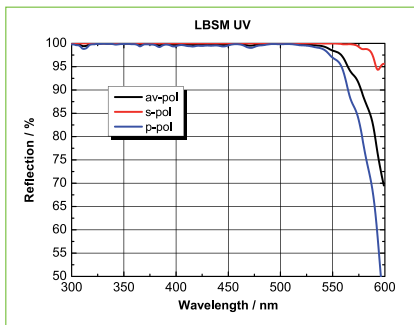
Dielectric-Coated Plane Mirrors LBSM-UV

- Dielectric coated plane mirror LBSM for laser applications
- Deflection mirror for 0-45° angles
- 1" (25.4 mm) diameter, 8 mm thickness
- Broadband reflection > 99 %
- Flatness λ/10
- Substrate: Borofloat
- Deflection mirror for 0-45° angles
- Reflection > 99 % for 300-550 nm for any polarization
- Damage threshold 500 kW / cm² (cw), 0.3 J/cm² (10 ns pulses)

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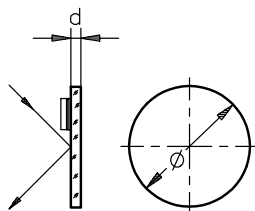
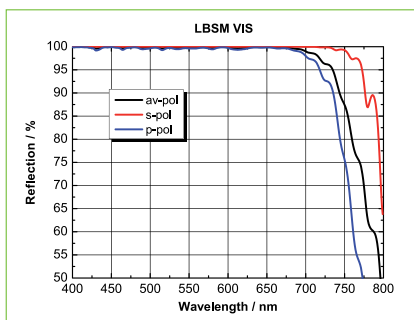
Dielectric-Coated Plane Mirrors LBSM-VIS

- Dielectric coated plane mirror LBSM for laser applications
- Deflection mirror for 0-45° angles
- 1" (25.4 mm) diameter, 8 mm thickness
- Broadband reflection > 99 %
- Flatness λ / 10
- Substrate: Borofloat
- Reflection > 99 % for 450-700 nm for any polarization
- Damage threshold 500 kW / cm² (cw), 1.0 J/cm² (10 ns pulses)

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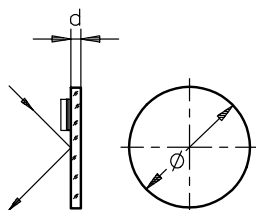
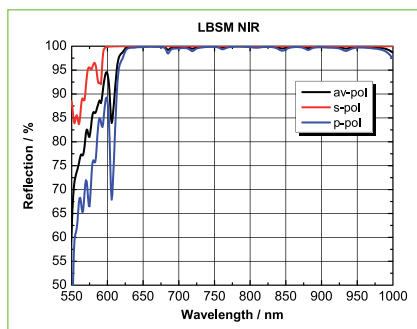
Dielectric-Coated Plane Mirrors LBSM-NIR

- Dielectric coated plane mirror LBSM for laser applications
- Deflection mirror for 45° angles
- 1" (25.4 mm) diameter, 8 mm thickness
- Broadband reflection > 99 %
- Flatness $\lambda/10$
- Substrate: Borofloat
- Reflection > 99 % for 700-900 nm
- Damage threshold 500 kW/cm² (cw), 10 J/cm² (28 ns pulses)

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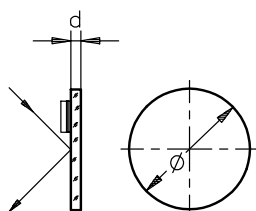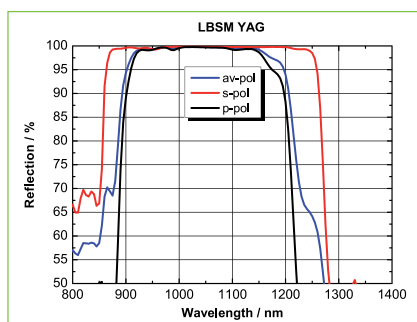
Dielectric-Coated Plane Mirrors LBSM-YAG

- Dielectric coated plane mirror LBSM for laser applications
- 1" (25.4 mm) diameter, 8 mm thickness
- Broadband reflection > 99 %
- Flatness $\lambda/10$
- Substrate: Borofloat
- Reflection > 99 % for 0-45° angles at 1064 nm
- Damage threshold 1 MW/cm² (cw), 5 J/cm² (10 ns pulses)

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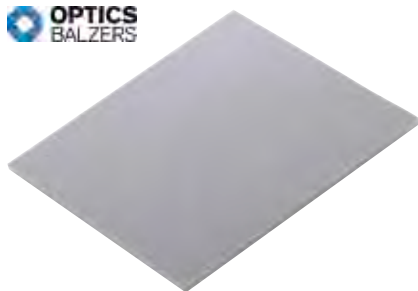Plane Mirror SILFLEX

- Mirrors for multimedia-, show- and projection systems
- No polarization and light effects
- Long-life
- High stability due to patented Sputter technique
- Substrate: heat-resistant Borosilicate glass
- SILFLEX-VIS: Extremely high reflection in the VIS ($R > 98\%$ for 420-680 nm, $\alpha = 0^\circ-45^\circ$)
- Special sizes up to 160 mm x 110 mm on request.

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
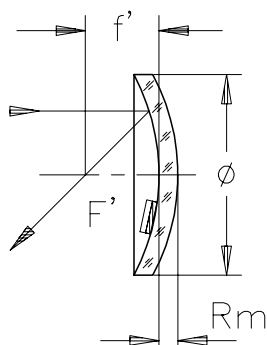
Front Surface Mirrors for Illumination Systems

- Designed for illumination systems
- Substrates are fabricated from low expansion material
- Aluminum coatings with a protective SiO_2 overcoating (RAL)
- Mirror radii = $2 \cdot f$
- The matching aspheric lenses are stated in the table.

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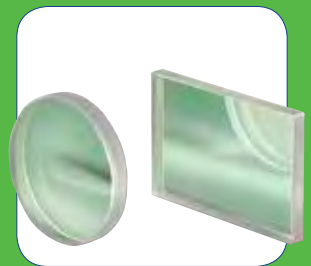Our quality criteria:

- *High-quality materials*
- *Continuous process monitoring*
- *Coatings for stringent demands in science and industry*
- *Laser-suitable coatings*
- *Expert consultation, from the development phase to the finished product*

LINOS Thin Film Coatings

From laser mirrors and filters to beamsplitters and anti-reflective coatings, the extensive range of Qioptiq coatings can meet your most stringent requirements.

If you can not find the coating that you need for your application among our standard coatings, please contact our experts, they will be happy to work with you in developing a special coating. With a wide range of production methods and many years of experience, our staff can offer a variety of solutions to find the best way to implement your designs. Our constant quality control ensures you reliability in planning, even for large series.

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Special Features:

To facilitate the development process we look forward to create preliminary designs, free of charge, based on your specifications.

Ideal Areas:

Laser optics; High-performance optics; Precision optics; Polarization optics; Special filters.

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Content

In our next Catalog this part (introduction; basics; information about measurement technology, environmental tests, job coating etc.) will also be available in English.

Price List for Standard Coatings 368

Anti-Reflective Coatings

Broadband Anti-Reflective Coating ARB 1 UV	369
Broadband Anti-Reflective Coating ARB 1	369
Broadband Anti-Reflective Coating ARB 1 IR	369
Broadband Anti-Reflective Coating ARB 2 UV	370
Broadband Anti-Reflective Coating ARB 2 VIS	370
Broadband Anti-Reflective Coating ARB 2 NIR	370
Broadband Anti-Reflective Coating ARB 2 IR	371
Broadband Anti-Reflective Coating ARB 3	371
Broadband Anti-Reflective Coating ARB 3 NIR	371
Broadband Anti-Reflective Coating ARB 4	372
Broadband Anti-Reflective Coating ARB 5	372
Broadband Narrowband Anti-Reflective Coating ARBS	373
High-Power Narrowband Anti-Reflective Coating ARHS UV	373
High-power Narrowband Anti-Reflective Coating ARHS	373
High-power doubleband anti-reflective coating ARHDS UV-VIS	374
High-power doubleband anti-reflective coating ARHDS VIS-NIR	374

Metallic Mirror Coatings

Reflective Aluminum Coating RAL-E UV	375
Reflective Aluminum Coating RAL VIS	376
Reflective Aluminum Coating RAL-E VIS	376
Reflective Aluminum Coating RAL-EE VIS	376
Front Surface Mirror Silver Coating RAGV	377
Rear Surface Mirror Silver Coating RAG	377
Reflective Gold Coating RAU	378
Reflective Gold Coating RAU-E	378

Dielectric Mirror Coatings

UV High-Power Laser Mirror Coating DLHS DUV	379
UV High-Power Laser Mirror Coating DLHS UV	379
High-Power Laser Mirror Coating DLHS VIS	379
High-power laser mirror coating DLHS Ti:Sa	380
IR High-Power Laser Mirror Coating DLHS NIR	380
Double Mirror Coating DLHD UV/VIS	380
Double Mirror Coating DLHD VIS/NIR	381
Broadband Mirror Coating DLB UV	381
Broadband Laser Mirror Coating DLB VIS	382
Broadband Mirror Coating DLB SBB	382

Beamsplitter Coatings

Partially Reflective Laser Mirror Coating TDLS	383
UV Broadband Beamsplitter Coating TBP UV	383
Broadband Beamsplitter Coating TBP VIS	384
Broadband Beamsplitter Coating TBP VIS NIR	384
NIR Broadband Beamsplitter Coating TBP NIR	384
IR Broadband Beamsplitter Coating TBP IR	385
Dielectric Broadband Beamsplitter Coating TBW UV	385
Dielectric Broadband Beamsplitter Coating TBW VIS	385
Broadband Hybrid Beam-splitter Coating TBW-H VIS	386
Polarizing Beamsplitter Cube Coating TSWP	386

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All specifications of our standard coatings can be found in our web-shop www.qioptiq-shop.com. Refer to the tab "Docs + Drawings" regarding the data sheets of the single coatings.

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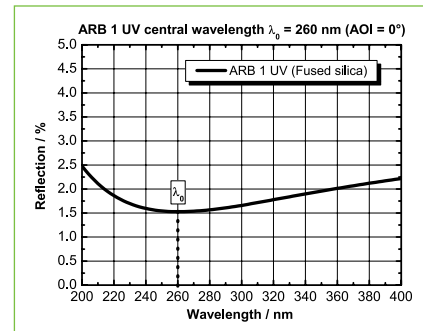
Broadband Anti-Reflective Coating ARB 1 UV

- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the UV spectral range*
- *Damage threshold: $H_{\infty} > 2 \text{ J/cm}^2$ at 308 nm with 10 ns laser pulses (s-on-1) at 10 Hz*
- *Design adaptable in the spectral range of $190 \text{ nm} \leq \lambda \leq 400 \text{ nm}$ and for larger angles of incidence*
- *Substrates: fused silica*

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
Broadband Anti-Reflective Coating ARB 1

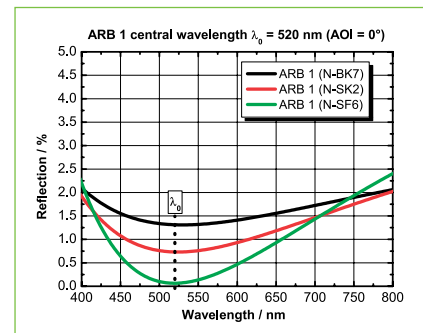- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission*
- *Residual reflection $0.3\% \leq R \leq 2\%$ depending on substrate for angles of incidence $0^\circ \leq \text{AOI} \leq 15^\circ$*
- *Damage threshold $H_{\infty} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *Design adaptable in the spectral range of $350 \text{ nm} \leq \lambda \leq 2400 \text{ nm}$ and for higher angles of incidence*
- *Substrates: N-BK7, fused silica or similar*

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
Broadband Anti-Reflective Coating ARB 1 IR

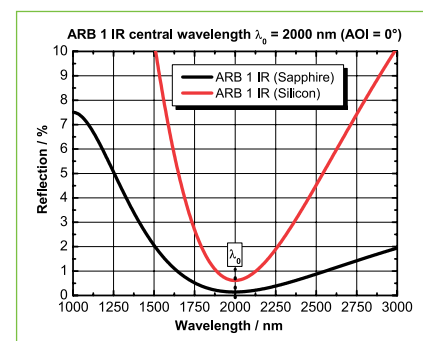- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the IR spectral range*
- *Damage threshold $H_{\infty} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz (Substrate: fused silica)*
- *Design adaptable in the spectral range of $1000 \text{ nm} \leq \lambda \leq 5000 \text{ nm}$ and for higher angles of incidence*
- *Substrates: sapphire, silicon*

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
Broadband Anti-Reflective Coating ARB 2 UV

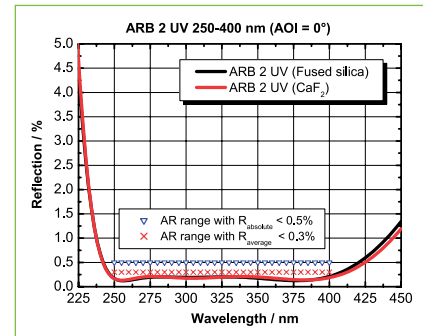- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the UV spectral range*
- *Damage threshold $H_{90} > 2 \text{ J/cm}^2$ at 266 nm with 9 ns laser pulses (s-on-1) at 10 Hz*
- *AR range, typical: 250-400 nm*
- *Design adaptable in the spectral range of $245 \text{ nm} \leq \lambda \leq 450 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, CaF_2*

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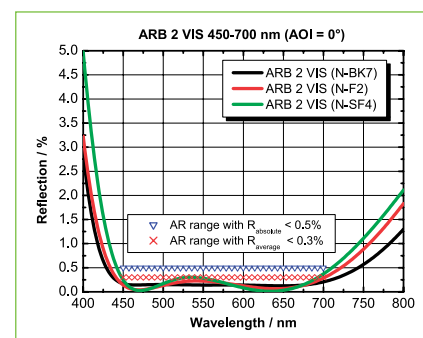
- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the visible spectral range*
- *Damage threshold $H_{90} > 10 \text{ J/cm}^2$ at 532 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *AR range typical: 450-700 nm*
- *Design adaptable in the spectral range of $350 \text{ nm} \leq \lambda \leq 950 \text{ nm}$ and for higher angles of incidence*
- *Substrates: N-BK7, fused silica or similar*

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
Broadband Anti-Reflective Coating ARB 2 NIR

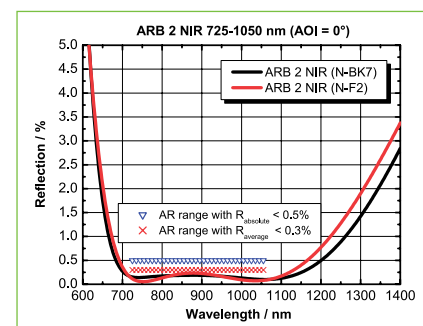- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the NIR spectral range*
- *Damage threshold: $H_{90} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *AR range typical: 725-1050 nm*
- *Design adaptable in the spectral range of $600 \text{ nm} \leq \lambda \leq 1400 \text{ nm}$ and for higher angles of incidence*
- *Substrates: N-BK7, fused silica or similar*

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
Broadband Anti-Reflective Coating ARB 2 IR

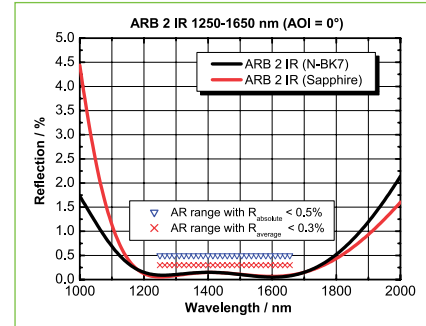- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the IR spectral range*
- *AR range typical: 1250-1650 nm*
- *Design adaptable in the spectral range of $1000 \text{ nm} \leq \lambda \leq 3000 \text{ nm}$ and for higher angles of incidence*
- *Substrates: sapphire, fused silica, N-BK7 or similar*

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
Broadband Anti-Reflective Coating ARB 3

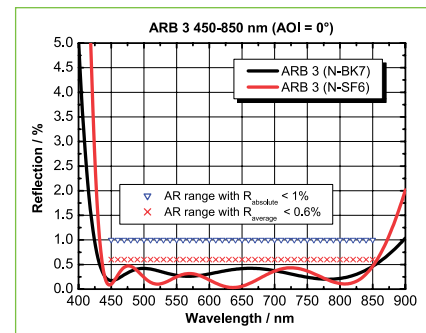- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission*
- *Damage threshold $H_{\infty} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *AR range typical: 450-850 nm*
- *Design adaptable in the spectral range of $260 \text{ nm} \leq \lambda \leq 950 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, N-BK7 or similar*

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
Broadband Anti-Reflective Coating ARB 3 NIR

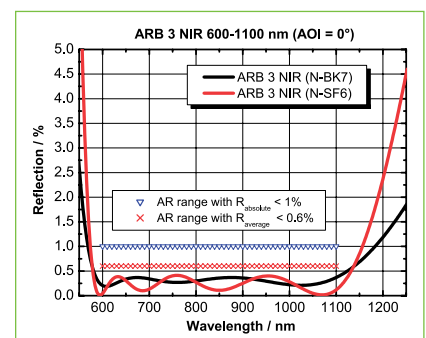- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the NIR spectral range*
- *Damage threshold: $H_{\infty} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *AR range typical: 600-1100 nm*
- *Design adaptable in the spectral range of $500 \text{ nm} \leq \lambda \leq 3000 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, N-BK7, high refractive glasses or similar*

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
Broadband Anti-Reflective Coating ARB 4

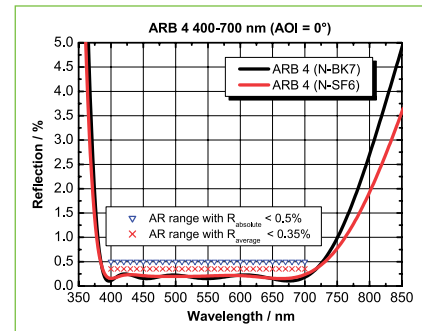- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission*
- *Damage threshold $H_{\infty} > 30 \text{ J/cm}^2$ for at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *AR range typical: 400-700 nm*
- *Design adaptable in the spectral range of $260 \text{ nm} \leq \lambda \leq 1500 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, N-BK7 or similar*

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
Broadband Anti-Reflective Coating ARB 5

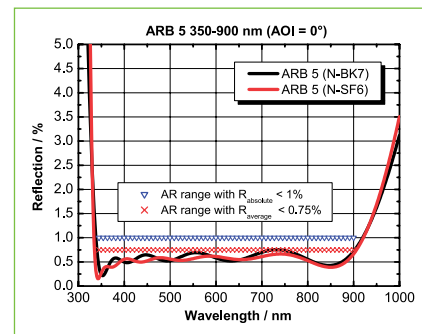- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission in an extremely broad spectral range*
- *Damage threshold $H_{\infty} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *AR range typical: 345-900 nm*
- *Design adaptable in the spectral range of $260 \text{ nm} \leq \lambda \leq 2200 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, N-BK7 or similar*

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You need special coatings for your optics? Please contact our experts for your individual request! To help you with precise facts please support us with the following data:

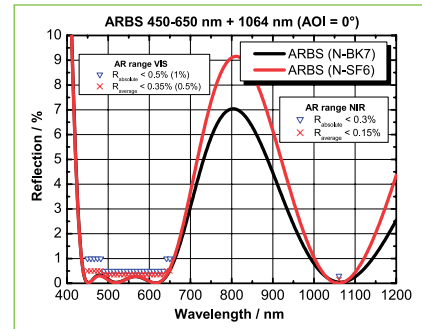
- **W**ellenlängenbereich
- **E**infallswinkel
- **S**ubstrat
- **P**olarisationseigenschaften
- **E**nergiedichte und weitere Umweltfaktoren
- *Wavelength range*
- *Angle of incidence*
- *Substrate*
- *Polarization specifications*
- *Energy and environmental facts*

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Broadband Narrowband Anti-Reflective Coating ARBS

- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for the visible spectral range in combination with a central wavelength in the NIR*
- *Damage threshold $H_{\infty} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *Design adaptable in the spectral range of $260 \text{ nm} \leq \lambda \leq 2200 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, N-BK7 or similar*
- *Please specify desired central wavelength λ_0 and wavelength range when ordering.*

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
High-Power Narrowband Anti-Reflective Coating ARHS UV

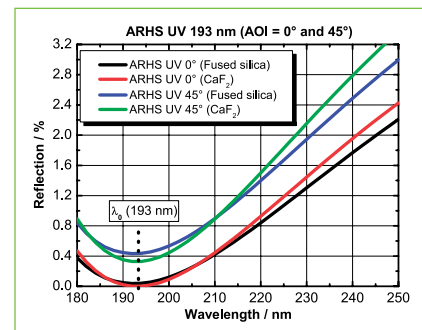- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission for high-power UV lasers*
- *Damage threshold $H_{\infty} > 1.8 \text{ J/cm}^2$ at 266 nm with 8.5 ns laser pulses (s-on-1) at 100 Hz*
- *Design adaptable in the spectral range of $180 \text{ nm} \leq \lambda \leq 400 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, CaF_2 or similar*
- *Please specify desired central wavelength λ_0 when ordering.*

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
High-power Narrowband Anti-Reflective Coating ARHS

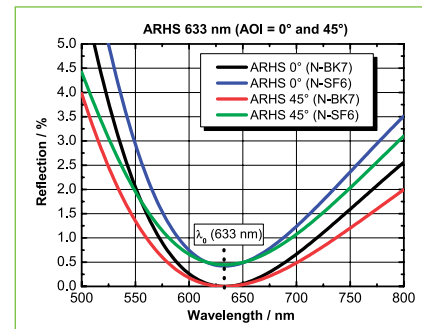- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission*
- *Design adaptable in the spectral range of $300 \text{ nm} \leq \lambda \leq 2200 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, N-BK7, high refractive glasses or similar*
- *Please specify desired central wavelength λ_0 as well as the AOI when ordering.*

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
High-power doubleband anti-reflective coating ARHDS UV-VIS

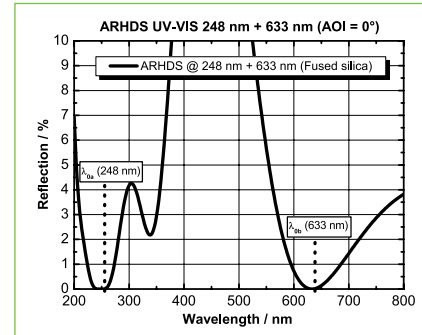- *Hard and scratch resistant anti-reflective coating to reduce reflection losses and increase transmission*
- *To use in the UV in combination with one alignment wavelength in the VIS*
- *Design adaptable in the spectral range of $180 \text{ nm} \leq \lambda \leq 950 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, CaF_2*
- *Please specify desired central wavelengths λ_0 when ordering.*

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
High-power doubleband anti-reflective coating ARHDS VIS-NIR

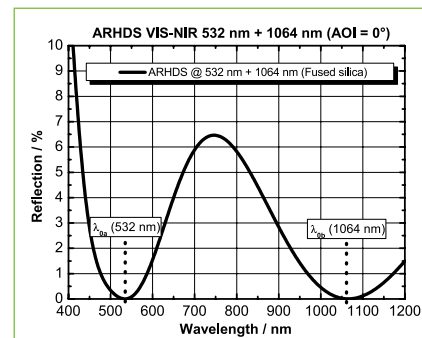- *Hard and scratch resistant anti-reflective coating for use in the NIR in combination with one alignment wavelength in the VIS to reduce reflection losses and increase transmission*
- *Design adaptable in the spectral range of $350 \text{ nm} \leq \lambda \leq 2200 \text{ nm}$ and for higher angles of incidence*
- *Substrates: fused silica, N-BK7, or similar*
- *Please specify desired central wavelengths λ_0 when ordering.*

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
Reflective Aluminum Coating RAL-E UV

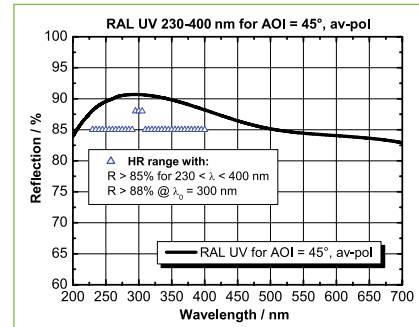- Standard aluminium reflective coating for mirrors for the UV spectral range
- High reflection within a broad spectral and angle-of-incidence range
- Guaranteed reflection $> 85\%$ in the range of $230\text{ nm} \leq \lambda \leq 400\text{ nm}$, $> 88\%$ at $\lambda_0 = 300\text{ nm}$ for $\text{AOI} = 45^\circ$, unpolarized
- Damage threshold $H_{90} > 200\text{ mJ/cm}^2$ at 193 nm with 30 ns laser pulses (s-on-1) at 10 Hz
- Design adaptable in the spectral range of $175\text{ nm} \leq \lambda \leq 450\text{ nm}$ and for higher angles of incidence
- Substrates: optical glasses, ceramic glasses, Zerodur, Borofloat, Pyrex, or similar

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Do you need special mirror coatings, anti-reflective coatings, beamsplitter coatings or polarization coatings for demanding applications?

We can provide expert consultation on all your requirements, from the development phase to production.

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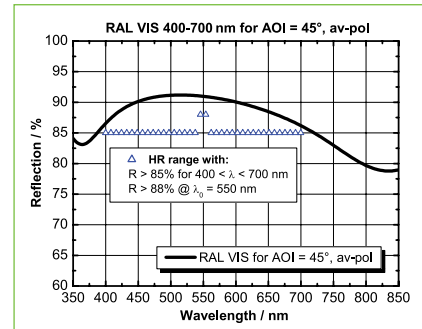
Reflective Aluminum Coating RAL VIS

- Standard aluminium reflective coating for mirrors in the VIS and IR spectral range up to 3 μm
- High reflection within a broad spectral and angle-of-incidence range
- Guaranteed reflection $> 85\%$ in the range of $400\text{ nm} \leq \lambda \leq 700\text{ nm}$, $> 88\%$ at $\lambda_0 = 550\text{ nm}$ for AOI = 45° , unpolarized
- Damage threshold $H_{\infty} > 1.5\text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Design adaptable in the spectral range of $350\text{ nm} \leq \lambda \leq 3000\text{ nm}$ and for higher angles of incidence
- Substrates: optical glasses, ceramic glasses, Zerodur, Borofloat, Pyrex, or similar

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
Reflective Aluminum Coating RAL-E VIS

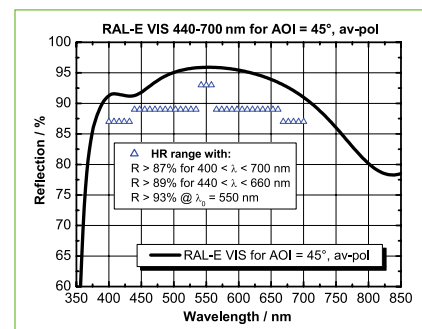- Enhanced reflective aluminium coating for mirrors for the VIS and NIR spectral range
- High reflection within a broad spectral and angle-of-incidence range
- Guaranteed reflection $> 89\%$ in the range of $400\text{ nm} \leq \lambda \leq 660\text{ nm}$, $> 93\%$ at $\lambda_0 = 550\text{ nm}$ for AOI = 45° , unpolarized
- Damage threshold $H_{\infty} > 1.5\text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Design adaptable in the spectral range of $450\text{ nm} \leq \lambda \leq 1000\text{ nm}$ and for other angles of incidence
- Substrates: optical glasses, ceramic glasses, Zerodur, Borofloat, Pyrex, or similar

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
Reflective Aluminum Coating RAL-EE VIS

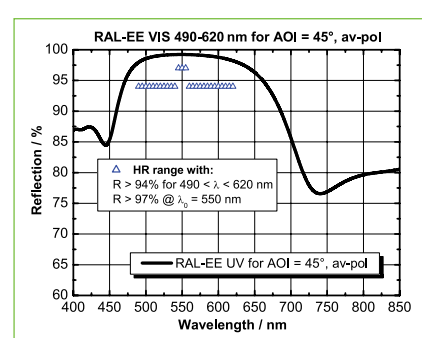- Enhanced high reflective aluminium coating for mirrors for the VIS and NIR spectral range
- High reflection within a broad spectral and angle-of-incidence range
- Guaranteed reflection $> 94\%$ in the range of $490\text{ nm} \leq \lambda \leq 620\text{ nm}$, $> 97\%$ at $\lambda_0 = 550\text{ nm}$ for AOI = 45° , unpolarized
- Damage threshold $H_{\infty} > 1.5\text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Design adaptable in the spectral range of $450\text{ nm} \leq \lambda \leq 1000\text{ nm}$ and for other angles of incidence
- Substrates: optical glasses, ceramic glasses, Zerodur, Borofloat, Pyrex, or similar

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
Front Surface Mirror Silver Coating RAGV

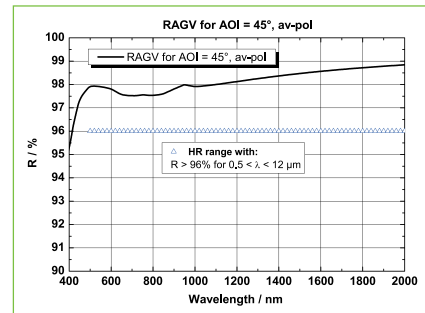- Front surface mirror silver coating for the VIS and IR spectral range up to 12 μm
- Constant high reflection within a broad spectral and angle-of-incidence range without absorption bands
- Guaranteed reflection ($0^\circ \leq \text{AOI} \leq 60^\circ$, unpolarized) $> 96\%$ in the range of $500 \text{ nm} \leq \lambda \leq 12 \mu\text{m}$
- Damage threshold $H_{10} > 1.5 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Substrates: optical glasses, ceramic glasses, Zerodur, Borofloat, Pyrex, or similar

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
Rear Surface Mirror Silver Coating RAG

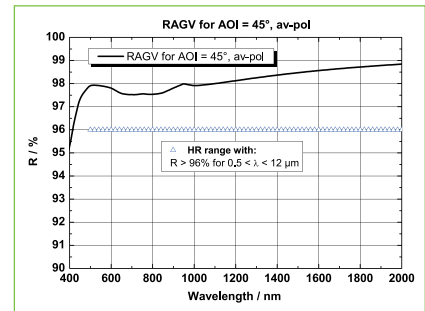- Metal film with the highest degree of reflection
- For rear surface mirrors
- Rear surface has a protective lacquer coating temperature stable from -40°C to $+100^\circ\text{C}$
- Guaranteed reflection $> 96\%$ in the range of $500 \text{ nm} \leq \lambda \leq 12 \mu\text{m}$
- Damage threshold $> 200 \text{ mJ/cm}^2$
- Substrates: optical glasses, ceramic glasses, or similar

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
Reflective Gold Coating RAU

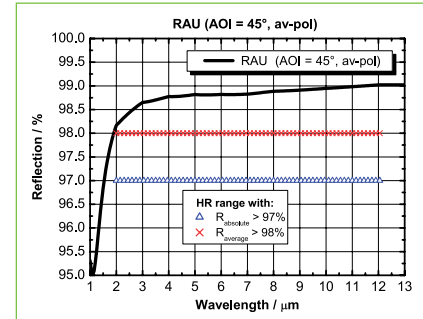- *Front surface mirror gold coating for the IR spectral range up to 20 μm*
- *Constant high reflection within a broad spectral and angle-of-incidence range without absorption bands*
- *Guaranteed reflection $> 97\%$ for $0^\circ \leq \text{AOI} \leq 60^\circ$ and $2 \mu\text{m} \leq \lambda \leq 20 \mu\text{m}$ (average polarization)*
- *Typical average residual reflection $> 98\%$ in the range of $2 \mu\text{m} \leq \lambda \leq 20 \mu\text{m}$ for $0^\circ \leq \text{AOI} \leq 60^\circ$, unpolarized*
- *Substrates: optical glasses, glass ceramic, diamond-turned Aluminium or Copper, or similar*

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
Reflective Gold Coating RAU-E

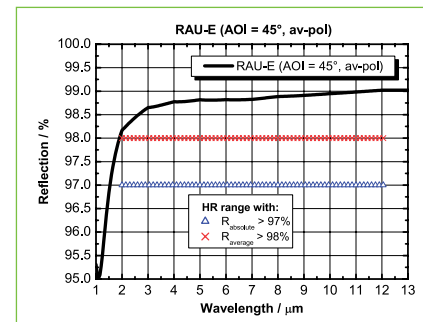- *Front surface mirror gold coating for the IR spectral range up to 25 μm with improved environmental characteristics*
- *Constant high reflection within a broad spectral and angle-of-incidence range without absorption bands*
- *Average reflection $> 98\%$ in the range of $0^\circ \leq \text{AOI} \leq 60^\circ$, unpolarized*
- *Abrasion resistant gold coating also on diamond turned or diamond milled aluminium*
- *Substrates: optical glasses, glass ceramic, diamond-turned Aluminium or Copper, or similar*

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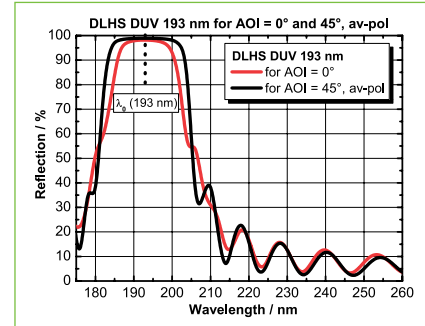
UV High-Power Laser Mirror Coating DLHS DUV

- Low-loss, hard and scratch resistant laser mirror coating for the DUV spectral range
- Guaranteed reflection > 96 % for central wavelength λ_0 for AOI = 0° or > 95 % for AOI = 45° (average polarization)
- Damage threshold $H_{10} > 5 \text{ J/cm}^2$ at 248 nm with 10 ns laser pulses (s-on-1) at 1 Hz
- Also available as dichroic
- Design adaptable in the spectral range of $180 \text{ nm} \leq \lambda \leq 260 \text{ nm}$ and for other angles of incidence
- Substrates: fused silica, CaF_2 , or similar
- Please specify desired central wavelength λ_0 and AOI (α) when ordering.


DLHS DUV for 193 nm (AOI = 0°) and DLHS DUV for 193 nm (AOI = 45°), unpolarized

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UV High-Power Laser Mirror Coating DLHS UV

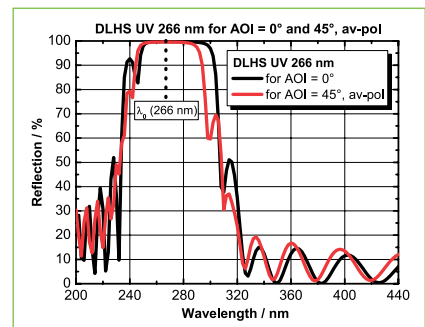- Low-loss, hard and scratch resistant laser mirror coating for the UV spectral range
- Damage thresholds $H_{10} > 5 \text{ J/cm}^2$ at 308 nm with 15 ns laser pulses (s-on-1) at 10 Hz
- Also available as dichroic
- Design adaptable in the spectral range of $240 \text{ nm} \leq \lambda \leq 450 \text{ nm}$ and for other angles of incidence
- Substrates: fused silica, CaF_2 or similar
- Please specify desired central wavelength λ_0 and AOI (α) when ordering.


DLHS UV for 266 nm (AOI = 0°) and DLHS UV for 266 nm (AOI = 45°), unpolarized

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High-Power Laser Mirror Coating DLHS VIS

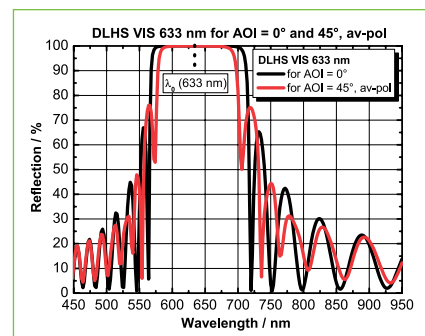- Low-loss, hard and scratch resistant laser mirror coating for the VIS spectral range
- Suitable for short pulse applications due to low dispersion of the coating ($\text{GDD} \sim 0 \text{ fs}^2$)
- Highest reflection within a broad spectral and angle-of-incidence range
- Guaranteed reflection > 99.8 % for central wavelength λ_0 for AOI = 0° or AOI = 45° (average polarization)
- Damage threshold $H_{10} > 10 \text{ J/cm}^2$ at 532 nm with 15 ns laser pulses (s-on-1) at 10 Hz
- Also available as dichroic
- Design adaptable in the spectral range of $350 \text{ nm} \leq \lambda \leq 950 \text{ nm}$ and for other angles of incidence
- Substrates: fused silica, N-BK7, or similar
- Please specify desired central wavelength λ_0 and AOI (α).

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
High-power laser mirror coating DLHS Ti:Sa

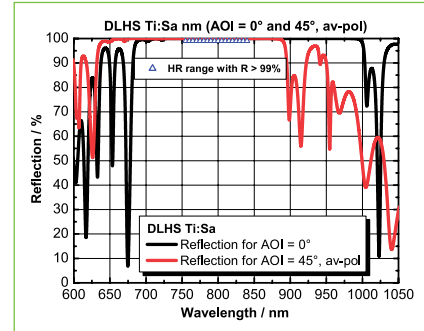- Low-loss, hard and scratch resistant laser mirror coating designed for Ti:Sapphire lasers
- Suitable for short pulse applications due to low dispersion of the coating (GDD ~ 0 fs²)
- Highest reflection within a broad spectral and angle-of-incidence range
- Guaranteed reflection > 99 % for the HR range for 0° ≤ AOI ≤ 45° (average polarization)
- Damage threshold $H_{\infty} > 10 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Substrates: fused silica and hard glass (e.g. Borofloat, Pyrex)


DLHS Ti:Sa for 755-840 nm (AOI = 0°) and DLHS Ti:Sa for 755-840 nm (AOI = 45°), unpolarized

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IR High-Power Laser Mirror Coating DLHS NIR

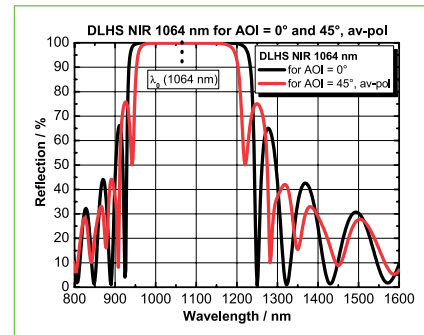- Low-loss, hard and scratch resistant laser mirror coating for the NIR spectral range
- Guaranteed reflection > 99.8 % for central wavelength λ_0 for AOI = 0° or AOI = 45° (average polarization)
- Damage threshold $H_{\infty} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Also available as dichroit
- Design adaptable in the spectral range of 750 nm ≤ λ ≤ 2200 nm and for other angles of incidence
- Substrates: fused silica, N-BK7, or similar
- Please specify desired central wavelength λ_0 and AOI (α) when ordering.

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
Double Mirror Coating DLHD UV/VIS

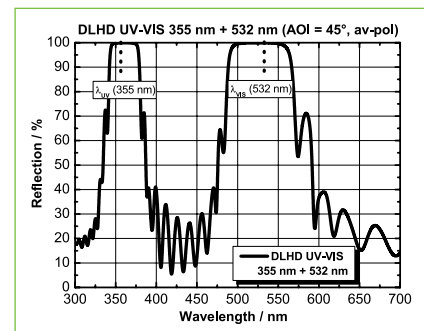- Low-loss, hard and scratch resistant laser mirror coating for two laser wavelengths, each in the UV and in the VIS spectral range
- Also suitable as beam steering mirror (AOI = 45°) or resonator end mirror (AOI = 0°)
- Guaranteed reflection > 99.8 % for laser wavelengths ($\lambda > 300 \text{ nm}$) in the UV and VIS spectral range for AOI = 0° or > 99.5 % for AOI = 45° (average polarization)
- Damage threshold $H_{\infty} > 30 \text{ J/cm}^2$ at 532 nm with 20 ns laser pulses (s-on-1)
- Design adaptable in the spectral range of 260 nm ≤ λ ≤ 1100 nm and for other angles of incidence
- Substrates: fused silica, N-BK7, or similar
- Please specify desired central wavelengths λ_1 and λ_2 and AOI (α) when ordering.

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
Double Mirror Coating DLHD VIS/NIR

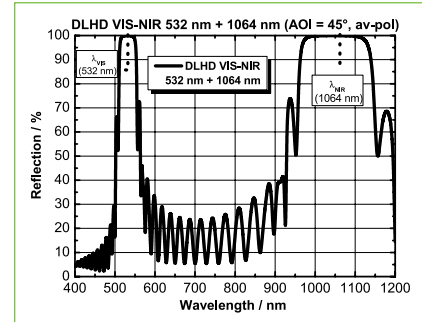- *Low-loss, hard and scratch resistant laser mirror coating for two laser wavelengths, each in the VIS and in the NIR spectral range*
- *Also suitable as beam steering mirror (AOI = 45°) or resonator end mirror (AOI = 0°)*
- *Guaranteed reflection > 99.8 % for laser wavelengths in the VIS and NIR spectral range or > 99.5 % for AOI = 45° (average polarization)*
- *Damage threshold $H_{\infty} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz*
- *Design adaptable in the spectral range of 350 nm $\leq \lambda \leq 1600$ nm and for other angles of incidence*
- *Substrates: fused silica, N-BK7, or similar*
- *Please specify desired central wavelengths λ_1 and λ_2 and AOI (α) when ordering.*

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
Broadband Mirror Coating DLB UV

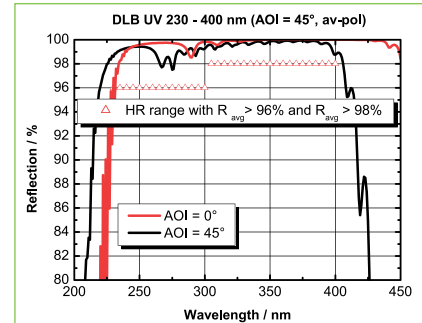- *Broadband and low-loss beam steering mirror coating for laser applications in the UV spectral range*
- *High reflection within a broad spectral and angle-of-incidence range*
- *$R_{\text{avg}} > 96\%$ at 230-300 nm ($0^\circ \leq \text{AOI} \leq 45^\circ$, unpolarized)*
- *$R_{\text{avg}} > 98\%$ at 300-400 nm ($0^\circ \leq \text{AOI} \leq 45^\circ$, unpolarized)*
- *Substrates: fused silica and hard glass (e.g. Borofloat, Pyrex)*

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
Broadband Laser Mirror Coating DLB VIS

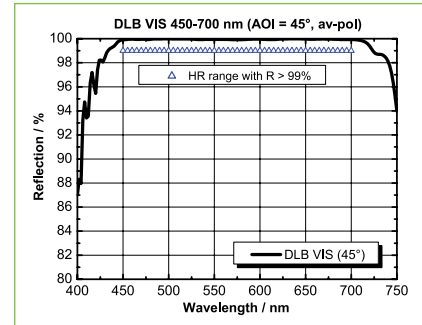- *Broadband and low-loss beam steering mirror coating for laser applications in the VIS spectral range*
- *Highest reflection within a broad spectral and angle-of-incidence range*
- *Guaranteed reflection > 99 % for the HR range for AOI = 45° (average polarization)*
- *Design adaptable in the spectral range of 390 nm ≤ λ ≤ 1600 nm and for other angles of incidence*
- *Substrates: fused silica and hard glass (e.g. Borofloat, Pyrex)*

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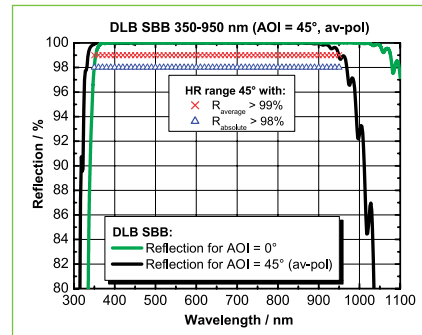
Broadband Mirror Coating DLB SBB

- *Extreme broadband and low-loss beam steering mirror coating for laser applications in the UV and NIR spectral range*
- *High reflection within a broad spectral and angle-of-incidence range*
- *Guaranteed reflection > 98 % for the HR range for AOI = 45° (average polarization)*
- *Design adaptable in the spectral range of 320 nm ≤ λ ≤ 1600 nm and for other angles of incidence*
- *Substrates: fused silica and hard glass (e.g. Borofloat, Pyrex)*

DLB SBB for 350-950 nm (AOI = 0°) and DLB SBB for 350-950 nm (AOI = 45°), unpolarized

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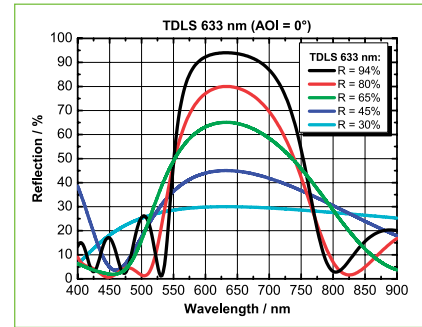
Partially Reflective Laser Mirror Coating TDLS

- *Low-loss, hard and scratch resistant laser mirror coating for laser applications in the UV, VIS and in the NIR spectral range*
- *Also suitable as output coupler (AOI = 0°) or as a plate beamsplitter*

- Any reflection ratio can be specified for the central wavelength λ_0

Damage thresholds:

- $H_{\infty} > 3 \text{ J/cm}^2$ at 248 nm with 20 ns laser pulses (s-on-1) at 10 Hz
 - $H_{\infty} > 5 \text{ J/cm}^2$ at 308 nm with 20 ns laser pulses (s-on-1) at 10 Hz
 - $H_{\infty} > 10 \text{ J/cm}^2$ at 532 nm with 15 ns laser pulses (s-on-1) at 10 Hz
 - $H_{\infty} > 30 \text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
 - Design adaptable in the spectral range of $180 \text{ nm} \leq \lambda \leq 2200 \text{ nm}$ and for other angles of incidence as well as for different polarization
 - Substrates: CaF_2 , N-BK7, fused silica, or similar
- Please specify desired central wavelength λ_0 and AOI (α) when ordering.

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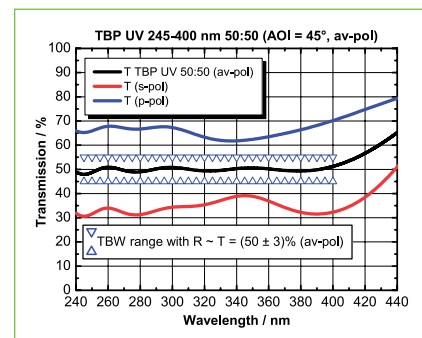
UV Broadband Beamsplitter Coating TBP UV

- *Low-loss, hard and scratch resistant beamsplitter coating for the ultra violet spectral range*
- *For splitting and combining of beam paths*

- Selectable transmission between $30 \% \leq T \leq 90 \%$
- Design adaptable within the spectral range from $240 \text{ nm} \leq \lambda \leq 500 \text{ nm}$, for other angles of incidence (AOI) and for certain polarizations
- Substrates: CaF_2 , fused silica, or similar

- Please specify desired wavelength range and angle of incidence when ordering.

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
Broadband Beamsplitter Coating TBP VIS

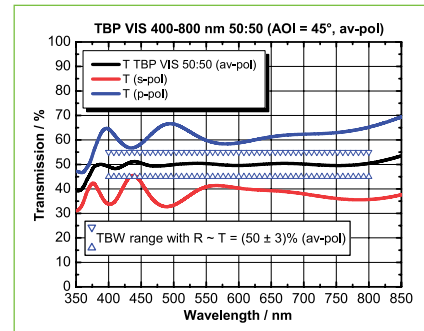- Low-loss, hard and scratch resistant beam-splitter coating for the visible spectral range
- For the splitting and combining of beam paths
- Selectable transmission between $30\% \leq T \leq 90\%$
- Design adaptable within the spectral range from $400\text{ nm} \leq \lambda \leq 1000\text{ nm}$, for other angles of incidence (AOI) and for certain polarizations
- Substrate: N-BK7, fused silica, or similar
- Please specify desired wavelength range and angle of incidence when ordering.

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
Broadband Beamsplitter Coating TBP VIS NIR

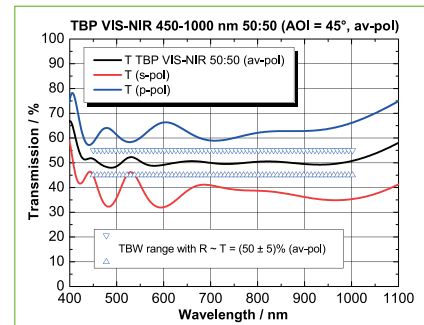- Low-loss, hard and scratch resistant beam-splitter coating for a broad spectral range from the VIS to NIR
- For splitting and combining of beam paths
- Selectable transmission between $30\% \leq T \leq 90\%$
- Design adaptable within the spectral range from $400\text{ nm} \leq \lambda \leq 1200\text{ nm}$, for other angles of incidence (AOI) and for certain polarizations
- Substrates: N-BK7, fused silica, or similar
- Please specify desired wavelength range and angle of incidence when ordering.

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
NIR Broadband Beamsplitter Coating TBP NIR

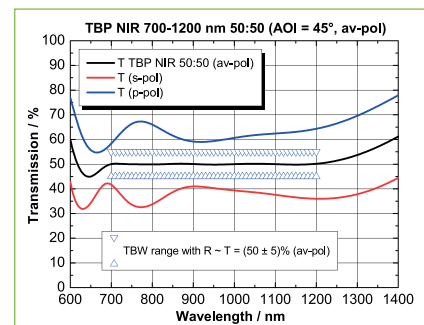- Low-loss, hard and scratch resistant beam-splitter coating for the near infrared spectral range
- For splitting and combining of beam paths
- Selectable transmission between $10\% \leq T \leq 80\%$
- Damage threshold $H_{10} > 10\text{ J/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Design adaptable within the spectral range from $600\text{ nm} \leq \lambda \leq 1400\text{ nm}$, for other angles of incidence (AOI) and for certain polarizations
- Substrates: N-BK7, fused silica, or similar
- Please specify desired wavelength range and angle of incidence when ordering.

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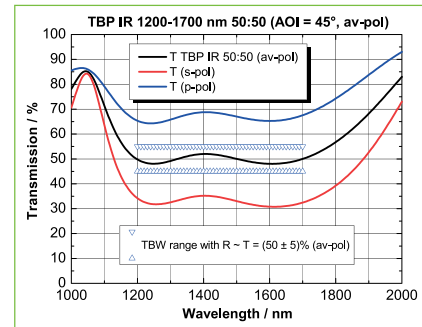
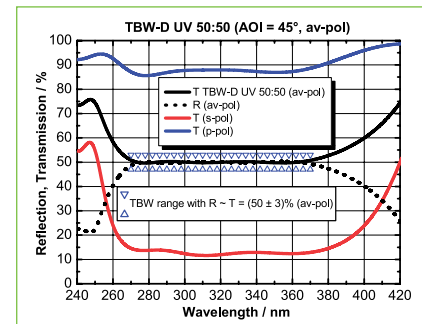
Dielectric Broadband Beamsplitter Coating TBW UV

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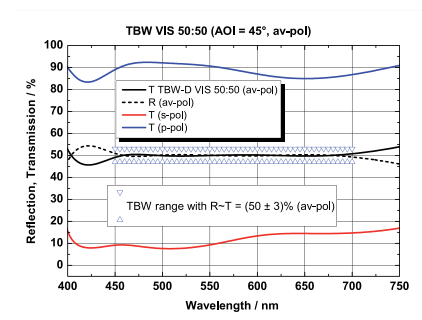
Dielectric Broadband Beamsplitter Coating TBW VIS

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
Broadband Hybrid Beam- splitter Coating TBW-H VIS

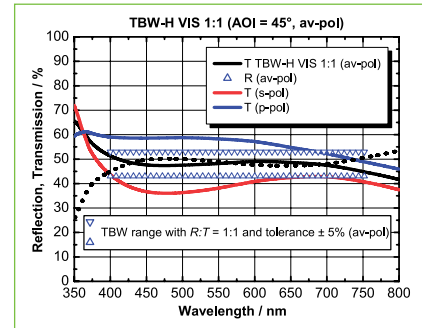- For beamsplitter cubes made from two cemented 90°-prisms
- For splitting and combining of beam paths
- Transmitted beam without beam displacement and beam deflection
- Applicable for low-power laser
- Selectable reflection between $30\% \leq T \leq 90\%$
- Damage threshold: limiting factor cemented system! Damage threshold $H_{\infty} > 100 \text{ mJ/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Design adaptable within the spectral range from $400 \text{ nm} \leq \lambda \leq 900 \text{ nm}$ and for certain polarizations
- Substrates: 90°-prisms from N-BK7, fused silica, or similar

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
Polarizing Beamsplitter Cube Coating TSWP

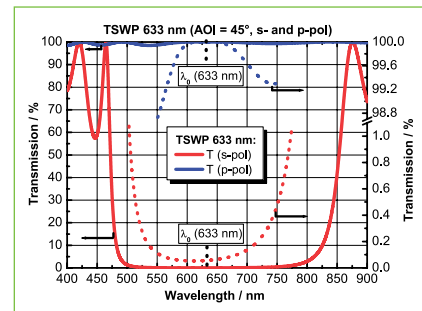- For polarizing beamsplitter cubes made from two cemented 90°-prisms
- Extreme splitting of polarization of the transmitted and reflected beam
- Guaranteed extinction ratio at central wavelength λ_0 for the transmitted beam $T_p : T_s > 1000 : 1$
- Damage threshold: limiting factor cemented system!
- Design adaptable within the spectral range from $400 \text{ nm} \leq \lambda \leq 2000 \text{ nm}$
- Substrates: glasses with a high refractive index like S-NBH 51, N-SF2, or similar

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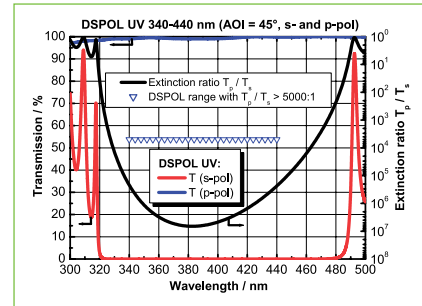
Thin Film Polarizer Coating DSPOL UV

- For polarizing cube beamsplitters made from two cemented 90°-triangle prisms for the ultra violet spectral range
- Extreme high splitting of polarization of the transmitted beam
- Guaranteed extinction ratio within the specified range of polarization (DSPOL-range) for the transmitted beam of $T_p : T_s > 5000 : 1$
- Damage threshold: limiting factor cemented system! Damage threshold $H_{\infty} > 20 \text{ mJ/cm}^2$ at 355 nm with 10 ns laser pulses (s-on-1)
- Design adaptable within the spectral range from $340 \text{ nm} \leq \lambda \leq 500 \text{ nm}$
- Substrates: N-BAK2 or similar

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
Thin Film Polarizer Coating DSPOL VIS

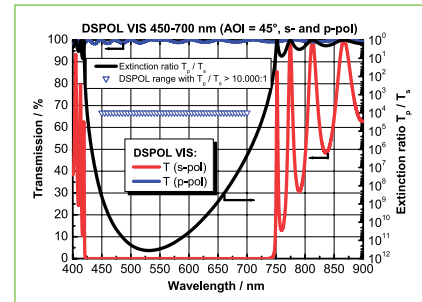- For polarizing cube beamsplitters made from two cemented 90°-prisms for the visible spectral range
- Extreme high splitting of polarization of the transmitted beam
- Guaranteed extinction ratio within the specified range of polarization (DSPOL-range) for the transmitted beam of $T_p : T_s > 10000 : 1$
- Damage threshold: limiting factor cemented system!
- Design adaptable within the spectral range from $400 \text{ nm} \leq \lambda \leq 850 \text{ nm}$
- Substrates: glasses with a high refraction index like S-NBH 51, N-SF4, or similar

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
Thin Film Polarizer Coating DSPOL VIS-NIR

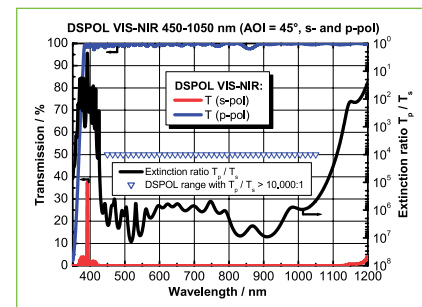- For polarizing cube beamsplitters made from two cemented 90°-prisms for the visible and the near infrared spectral range
- Extreme high splitting of polarization of the transmitted beam within a broad spectral range from the visible up to the near infrared range
- Guaranteed extinction ratio within the specified range of polarization (DSPOL-range) for the transmitted beam of $T_p : T_s > 10000 : 1$
- Damage threshold: limiting factor cemented system! Damage threshold $H_{\infty} > 100 \text{ mJ/cm}^2$ at 1064 nm with 11 ns laser pulses (s-on-1) at 10 Hz
- Design adaptable within the spectral range from $400 \text{ nm} \leq \lambda \leq 1200 \text{ nm}$
- Substrates: S-NBH 51, N-SF4, or similar

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
Thin Film Polarizer Coating DSPOL IR

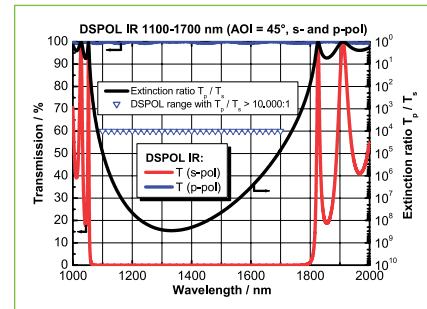- For polarizing cube beamsplitters made from two cemented 90°-prisms for the infrared spectral range
- Extreme high splitting of polarization of the transmitted beam
- Guaranteed extinction ratio within the specified range of polarization (DSPOL-range) for the transmitted beam of $T_p : T_s > 10000 : 1$
- Damage threshold: limiting factor cemented system!
- Design adaptable within the spectral range from $750 \text{ nm} \leq \lambda \leq 2200 \text{ nm}$
- Substrates: S-NBH 51, N-SF4, or similar

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Our quality criteria:

- *High-quality materials*
- *High-quality polarization coatings*
- *Low wave front distortion*
- *The narrowest possible production tolerances*

Polarization Optics

Qioptiq offers an extensive line of polarizers, polarizing filters and wave plates made of high-quality materials for the best possible transmission.

Complex coatings ensure high extinction ratios and broadband solutions. Our narrow optical production tolerances ensure minimal wave front distortion.

And of course, our polarization optics are available in Microbench compatible mounts.

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Special Features:

Optionally mounted for LINOS bench systems

Ideal Areas:

Polarization rotation; voltage measurement; polarization-dependent beam splitting.

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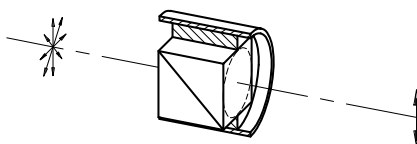
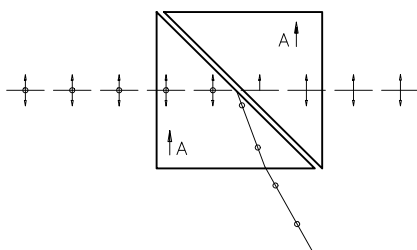
Maximum angular polarized field as a function of wavelength:

- a) full angular field
b) symmetrical half-angular field

	214 nm	298 nm	361 nm	404 nm	589 nm	1041 nm
a)	12.7°	9.9°	8.8°	8.5°	8.0°	7.5°
b)	1.7°	3.8°	4.3°	3.8°	2.7°	2.0°

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Glan-Taylor Prism Polarizers

- Polarizing cube composed of two air-spaced calcite prisms with their optical axes aligned mutually parallel
- Extremely large extinction ratios
- Virtually no deviation of extraordinary ray
- Mounted, outer diameter fits to Microbench

- Uncoated
- Length / aperture ratio = 0.85
- Surface flatness : $\lambda/4$ @ 633 nm over the clear aperture
- Surface quality : 40-20 scratch dig, i.e. 5/2x0.16 (DIN ISO 10110)
- Transmission range 250-2300 nm, transmission (uncoated):
25-30 % at 250 nm
40-45 % at 300 nm
65-70 % at 400 nm
85-87 % beyond 500 nm

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
These prism polarizers are also available without mount on request. Please write to sales.gioptiq@excelitas.com or give us a call +49(0).551 6935-0. We are pleased to perform your requirements.

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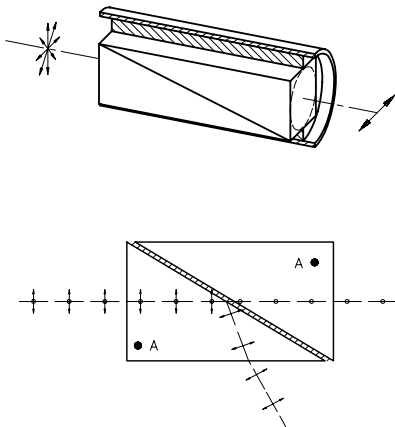Glan-Thompson Prism Polarizers


- Composed of two cemented calcite prisms with their optical axes aligned parallel to their hypotenuse faces
- Extremely large extinction ratios
- Angular fields are larger than those of Glan-Taylor prisms
- Angular field symmetric about prism assembly axis
- Virtually no deviation of extraordinary ray
- Mounted, fits to Microbench
- Uncoated
- Damage thresholds: CW lasers: 1 W/cm²; Pulsed lasers: 5 W/cm²
- Useable angular polarized field symmetrical with prism axis 14° - 16°
- Length/aperture ratio = 2.5
- Surface flatness : $\lambda/4$ @ 633 nm over the clear aperture
- Surface quality : 40-20 scratch dig, i.e. 5/ 2x0.16 (DIN ISO 10110)
- Transmission range 350-2300 nm, transmission (uncoated):
35-40 % at 350 nm
65-70 % at 400 nm
80-88 % beyond 500 nm
- Optionally available unmounted.

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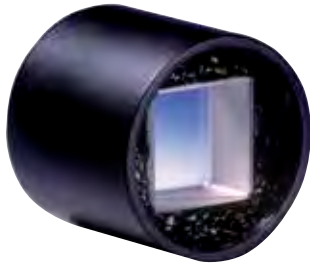¹⁾ Material classifications: S: striation free, i.e., no detectable variations in refractive index; no striations discernible under schlieren testing; A: slight variations in refractive index, i.e., exhibiting minor striations, just barely discernible under schlieren testing

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


Rochon Prism Polarizer

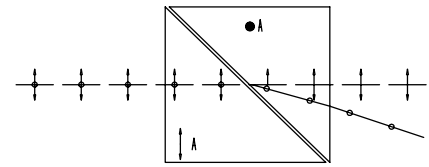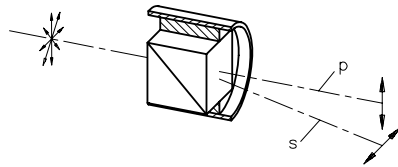- Consists of two calcite prisms
- Extremely large extinction ratios
- Ordinary ray is achromatic
- Mounted, fits to Microbench
- Uncoated
- Maximum deviation of the ordinary ray is better than 3'
- Parallelism of end surfaces $\lambda/4$
- Surface flatness : $\lambda/4$ @ 633 nm over the clear aperture
- Surface quality : 40-20 scratch dig, i.e. 5/ 2x0.16 (DIN ISO 10110)
- Transmission (uncoated): > 85 % at 350-2500 nm
- Optionally unmounted available.

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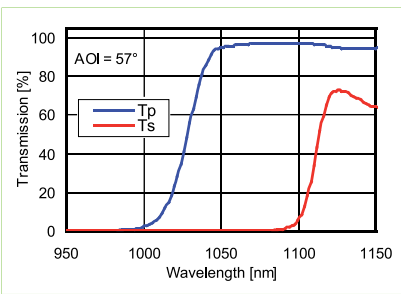
Fused Silica Brewster Polarizer


- For complete polarization of light
- Rectangular design
- Fused silica
- Extinction ratio > 500:1 at 1064 nm
- Angel of incidence = 57° ± 2°
- Damage thresholds (1064 nm): > 30 J/cm² at 10 ns; > 120 J/cm² at 160 ns; > 1 MW/cm² at cw

Brewster-Polarisator aus Quarzglas Brewster Polarizer fabricated from fused silica

Größe lxb (mm)	Dicke d (mm)	Toleranz lxb (mm)	Toleranz d (mm)	Planität beidseitig (λ)	Parallelität (mm)	Sauberkeit nach ISO 10110	Part No.
Dimensions lxb (mm)	Thickness d (mm)	Tolerance lxb (mm)	Tolerance d (mm)	Flatness Both Surfaces (λ)	Parallelism (mm)	Surface Quality ISO 10110	Part No.
14.3x28.6	3.2	±0.1	±0.1	1/10	20"	5x0.025	G390232000

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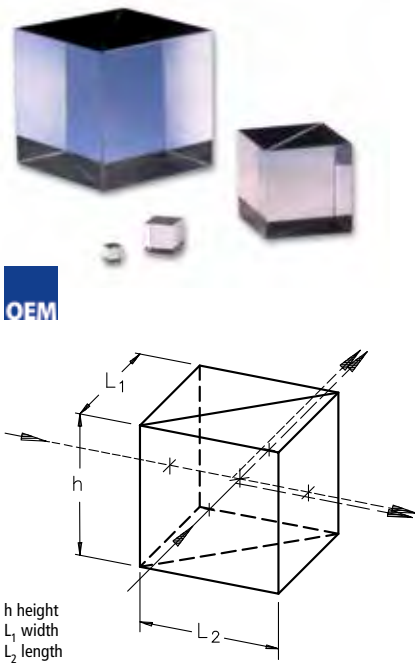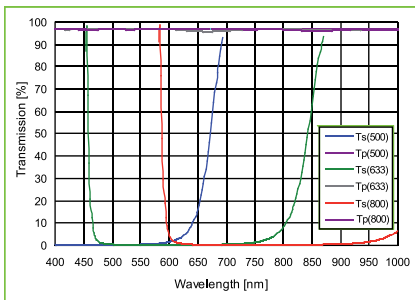Polarizing Beamsplitter Cubes

- For 450-550 nm, 550-700 nm, and 700-900 nm (TSWP 800)
- Over-all transmission of the cubes $T_p \geq 90\%$, typical $T_p \geq 95\%$
- Broadband ARB2 coatings
- 8' angular beam deflection tolerance
- Damage thresholds: > 2 kW/cm² cw (488/514 nm)
- Damage thresholds: > 200 mJ/cm² for pulses of 10 ns (1064 nm)
- Damage thresholds: > 28 mJ/cm² for pulses of 280 fs (850 nm, 20 Hz)

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
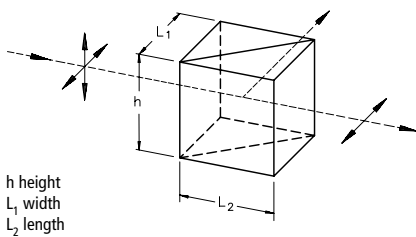
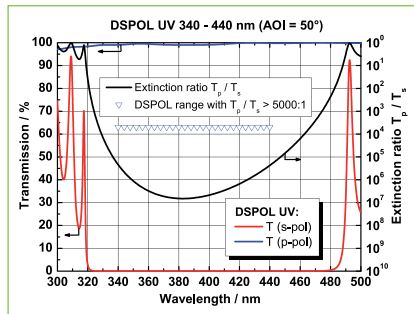
Thin Film Polarizers

- *Special beamsplitter coatings yield highly polarized transmitted beams*
- *No angular or lateral deflection of transmitted beams*
- *Fits to Microbench with clamp holder*
- *Broadband coated with ARB2*
- *Combine extreme spectral bandwidth with superb extinction ratio and transmittance*


- $T_p > 90\%$
- *Extinction ratio transmission $> 10,000:1$ ($\Phi_{Tp} : \Phi_{Ts}$) at $\pm 2^\circ$*
- *Less than $\lambda/4$ wavefront distortion at 633 nm*
- *8' angular beam deflection tolerance*
- *Damage thresholds: $> 2 \text{ kW/cm}^2 \text{ cw}$ (488/514 nm)*
- *Damage thresholds: $> 200 \text{ mJ/cm}^2$ for pulses of 10 ns (1064 nm)*
- *Damage thresholds: $> 28 \text{ mJ/cm}^2$ for pulses of 280 fs (850 nm, 20 Hz)*

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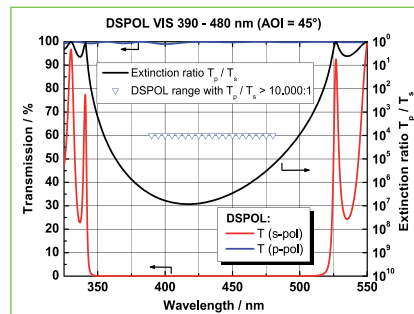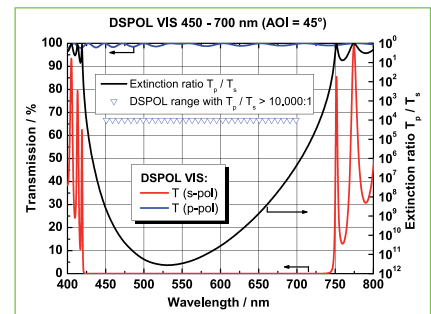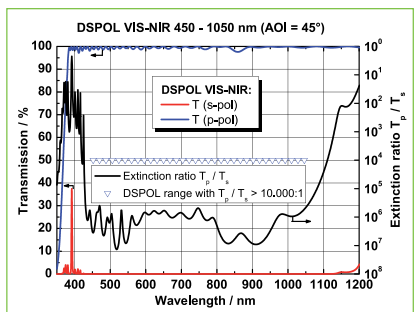
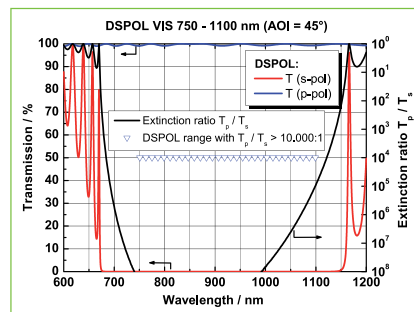
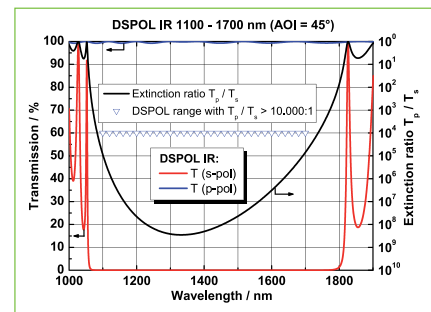
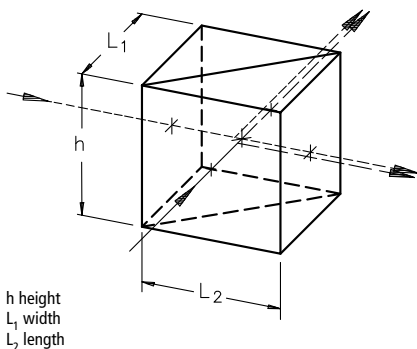
Optical Contacted Polarizing Beamsplitter Cubes

These polarizing beamsplitter cubes are derived from a unique combination of ultraprecise fabrication capabilities and epoxy-free bonding technology which results in a zero-bondline thickness.

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

- High-energy laser-line and broadband designs
- Optical bonding for epoxy-free optical paths
- Convenient 90° beam separation

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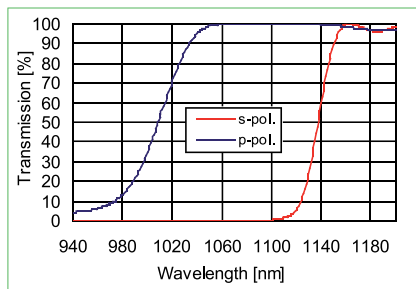
- Surface Flatness $\lambda/10$
- Transmitted Beam Deviation < 0.9 mrad
- Average transmission > 95 %
- Fabricated from fused silica

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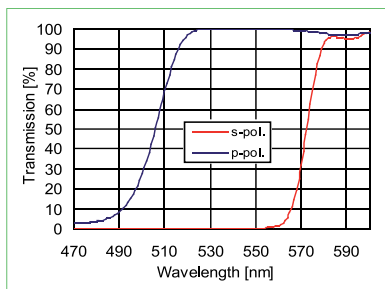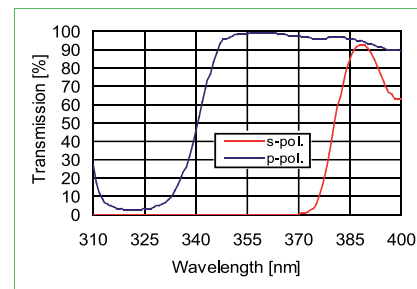Take advantage of more than 130 years of experience in the manufacture of precision optics and precision mechanical components and systems.

Früher
In former times

Heute
Today

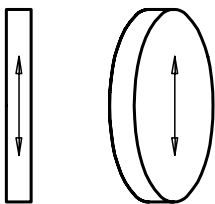
Montage und Mechanikfertigung in
Göttingen um 1910
*Mounting and mechanics production in
Goettingen about 1910*


Einstellung der Parameter an einer unserer
Coatinganlagen
*Setting of parameters at one of our
coating units*

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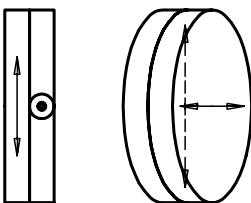Retardation Plates

- Made of crystal quartz
- Antireflective coating ($R < 0.2\%$ at central wavelength)
- Wavefront distortion: $\lambda/8$ (at 632.8 nm)
- Surface quality: 2×0.063 (20/10 scratch and dig)
- Parallelism better than $1''$
- Retardation tolerance $\lambda/500$ (typical)

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
Type S retarders are among the higher-order retarders, because the overall thickness of the plate is not just $\lambda/2$ or $\lambda/4$, but also a whole-number multiple of λ . This makes the plates easier to manufacture and handle.

With Type D, this is compensated through combination with a second retarder with its axes rotated in relation to those of the first plate. The resulting retardation is thus no more than the desired fraction of λ (without the whole-number multiple). The advantage of Type D retarders lies in their low temperature sensitivity.

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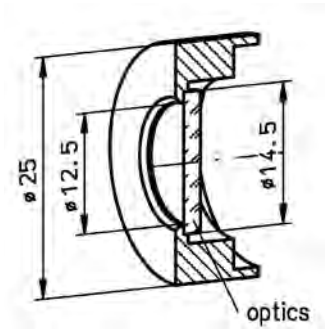
Achromatic Retarders

Practical and versatile: In contrast to simple wave plates the combination of different crystal plates of magnesium fluoride and quartz-defined enables defined delays for broad wavelength ranges. Therefore one component can be used for several different laser lines.


- Retardations: $\lambda/2$ and $\lambda/4$
- Three different broad wavelength ranges
- Mounted versions fit to Microbench
- Broadband
- Cemented
- Transmission > 86 %
- Uncoated; coating available on request
- Surface flatness: $\lambda/4$ at 633 nm (over clear aperture)
- Surface quality: 40-20 (5/ 2x0.16 according to DIN ISO 10110)
- Wavefront distortion: $\lambda/2$ up to 1λ
- Damage threshold: > 2 kW/cm² cw (488 /514 nm)
- Damage threshold: > 200 mJ/cm² for pulses of 10 ns (1064 nm)

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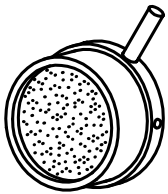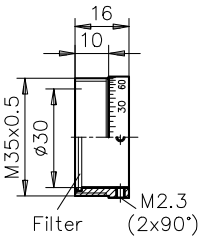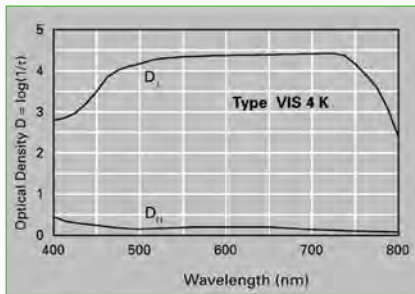
Polarization Filter Type VIS 4 K

- Linearly polarizes dichroic sheet polarizer
 - Ideally suited for demonstrations or testing, stress analysis, and for continuously attenuating linearly polarized light at constant clear aperture
 - Strikes a good compromise between high transmittance and high extinction ratio
 - Unmounted and mounted (two types of mount) available
- Transmission for unpolarized white light ≈ 32 %
 - Resultant transmittance of two filters, E-vector direction parallel (DP) > 20 %
 - Resultant transmittance of two filters, E-vector directions crossed (DS) < 0.005 %
 - Extinction ratio 1 : 4,000
 - Degree of linear polarization > 99.97 %
 - Max. long-term temperature ≈ 70 °C

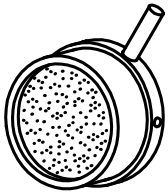
Mounted types in:

- Rotary insert with scale 0° to 180° for Microbench mounting plates G06 1045 000 and G06 1024 000
- or Mount Ø 25 mm with adjusting lever for Microbench components with inner Ø 25 mm.

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
Qioptiq uses glass from established manufacturers in the production of precision optics components. This ensures the consistent high quality of our polarization optics.

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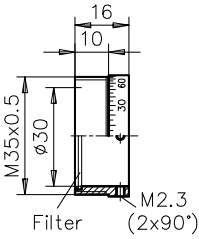Retardation Sheets $\lambda/4$

Retardation sheets are achromatic sheets of 0.3 mm thickness. These sheets work according to conventional Gypsum- or Mica plates. Quarter-waveplate sheet especially suitable in the wavelength range of 480-640 nm. Please combine with a linear polarizer (in an angle of 45°) to generate circular polarized light.

- Unmounted and mounted (two types of mount) available

- Mounted types in:
- Rotary insert with scale 0° to 180°, 5° scaling for Microbench mounting plates G06 1045 000 and G06 1024 000
 - or Mount Ø 25 mm with adjusting lever for Microbench components with inner Ø 25 mm.

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Polarization Filter 80 in Rotary Mount

Thin sheet polarizer without plastic protective film, directly cemented between plano cover glass

- Ideal for stress analysis and polarimetry
- Sizable free aperture
- Linear polarizing
- Mount with 1° scaling
- Mean transmittance of single filter for unpolarized white light $\approx 42\%$
- Resultant transmittance of two filters, E-vector direction parallel (D_p) $> 38.5\%$
- Resultant transmittance of two filters, E-vector directions crossed (D_d) $< 0.01\%$
- Extinction ratio 1:3,500
- Degree of linear polarization $> 99.5\%$
- Max. long-term temperature $\approx 70^\circ\text{C}$
- Rotary mount with 1°-scale
- 2.6 mm thick
- Free aperture: 78 mm
- Mounting options: central M6 tapped hole, 2xM4 / 50 mm centers

Polfilter 80 in Drehfassung Polarization filter 80 in rotary mount

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Polarization filter 80 in rotary mount	G036325000

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Our quality criteria:

- *Sophisticated design*
- *Many years of experience*
- *Latest optical, mechanical and mount technology*

Laseroptics

Benefit from our many years of experience in the development of optical systems for laser material processing.

Our broad selection of beam expansion systems and F-Theta-Ronar lenses meets even the most stringent demands. This comprehensive range covers everything from fixed beam expanders to variable and motorized beam expanders. The high numerical aperture of LINOS laser lenses has gained them wide popularity.

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Special Features:

F-Theta Ronar objectives: optical glass, fused silica and telecentric lenses available.

Ideal Areas:

Marking, structuring, perforating, drilling and more for laser material processing.

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Beam Expanders

<i>LINOS Variable Magnification Beam Expanders</i>	416
<i>LINOS Motorized Beam Expanders</i>	418
<i>Laser Beam Expander System 4x and 7x</i>	420
<i>Laser Beam Expander System 10x</i>	420

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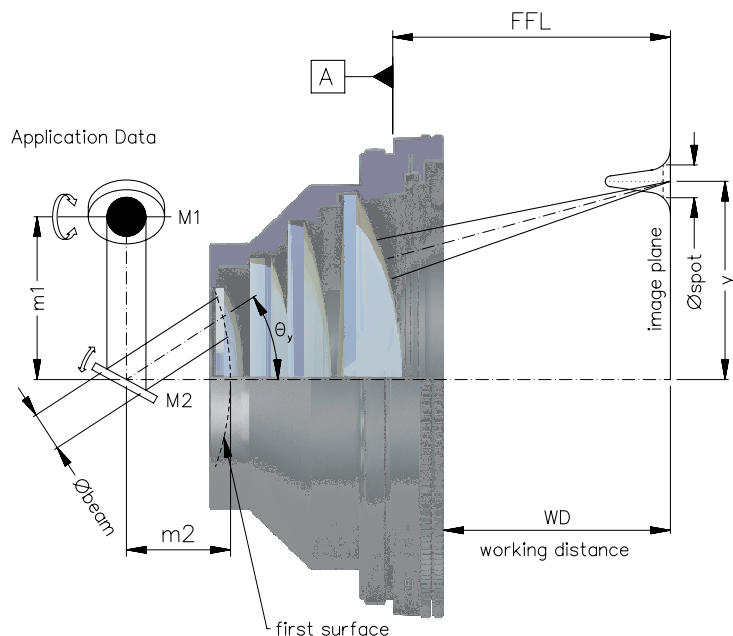
LINOS F-Theta-Ronar Lenses 340-360 nm

- Fused-silica designs
 - Telecentric versions available
 - Focal lengths ranging from 70 mm to 255 mm, tolerance $\pm 1\%$
 - Screw thread M85x1
- Transmission $\geq 96\%$ with good performance in VIS-range
 - Laser-damage threshold coating up to 4 J/cm^2 at 355 nm, 6 ns, 100 Hz
 - Includes interchangeable fused-silica protective glasses

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All LINOS F-Theta-Ronar lenses can be used with enlarged beam diameters and different mirror distances. Accordingly the scan fields and spot size diameters will be changed. Please feel free to send us your request.

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LINOS F-Theta-Ronar Lenses 515-540/532 nm

- Fused-silica and optical-glass designs
- Telecentric versions available
- Focal lengths ranging from 100 mm to 420 mm, tolerance $\pm 1\%$
- Screw thread M85x1
- Transmission $\geq 96\%$ with good performance in VIS-range
- Laser damage threshold coating up to 20 J/cm² at 532 nm, 6 ns and up to 0.7 J/cm² at 515 nm, 285 fs
- Includes interchangeable protective glasses

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Fused silica lenses

Qioptiq has developed a range of sophisticated F-Theta-Ronar scan lenses made of fused silica for high-power and short-pulse laser material processing. Fused-silica lenses provide minimized thermal focus shift and higher resistance when working at high power density. A specially developed coating achieves very low reflection and absorption values.

LINOS F-Theta-Ronar 515-540/532 nm, Quarzglas LINOS F-Theta-Ronar 515-540/532 nm, Fused Silica

Nominal-brennweite (mm)	Max. Scanfeld (mm ²)	Scanwinkel $\pm \Theta_{x,y}$ max. (°)	Strahlenbündel-Durchmesser begrenzt bei 1/e ² \varnothing_{beam} (mm)	Bildpunkt-Durchmesser 1/e ² \varnothing_{spot} (µm)	Spiegelabstände (m1/m2) (mm)	Arbeitsabstand (WD) (mm)	Schutzglas	Artikel-Nr.
Nominal focal length (mm)	Maximum scan field (mm ²)	Scan angle $\pm \theta_{x,y}$ max. (°)	Beam diameter truncated at 1/e ² \varnothing_{beam} (mm)	Spot diameter at 1/e ² \varnothing_{spot} (µm)	Mirror distances m1/m2 (mm)	Working distance WD (mm)	Protective glass	Part No.
100 telecentric	43 x 43	± 12.2	14	9	17/28	138.5	PG13	4401-547-000-21
167 telecentric	86 x 86	± 15.4	14	12	17/33	215.5	PG21	4401-517-000-21
255	170 x 170	± 19.3	10	25	13/30	318.1	PG13	4401-496-000-21

LINOS F-Theta-Ronar 515-540/532 nm, Optisches Glas LINOS F-Theta-Ronar 515-540/532 nm, Optical glass

Nominal-brennweite (mm)	Max. Scanfeld (mm ²)	Scanwinkel $\pm \Theta_{x,y}$ max. (°)	Strahlenbündel-Durchmesser begrenzt bei 1/e ² \varnothing_{beam} (mm)	Bildpunkt-Durchmesser 1/e ² \varnothing_{spot} (µm)	Spiegelabstände (m1/m2) (mm)	Arbeitsabstand (WD) (mm)	Schutzglas	Artikel-Nr.
Nominal focal length (mm)	Maximum scan field (mm ²)	Scan angle $\pm \theta_{x,y}$ max. (°)	Beam diameter truncated at 1/e ² \varnothing_{beam} (mm)	Spot diameter at 1/e ² \varnothing_{spot} (µm)	Mirror distances m1/m2 (mm)	Working distance WD (mm)	Protective glass	Part No.
100	58 x 58	± 17.7	6	15	16/12	90.8	PG8	4401-304-000-21
100 telecentric	54 x 54	± 15.6	15	7	20/32	126.6	PG7	4401-461-000-21
160	98 x 98	± 17.7	10	16	16/12	176.1	PG8	4401-305-000-21
250	154 x 154	± 17.7	20	12	22/24	288.5	PG7	4401-289-000-20
330	204 x 204	± 17.7	14	23	18/24	389	PG7	4401-485-000-21
420	275 x 275	± 18.7	15	27	17/28	494.6	PG7	4401-489-000-21

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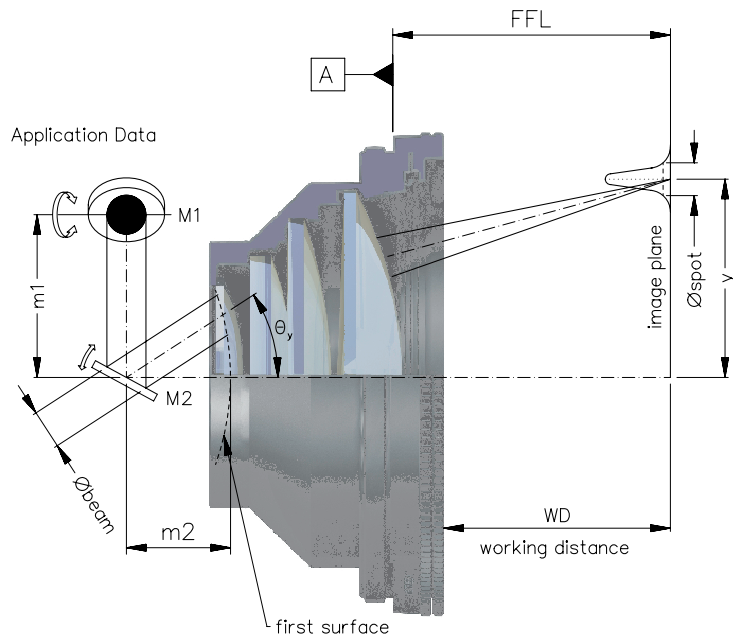
LINOS F-Theta-Ronar Lenses 940-980 nm

- Focal lengths ranging from 100 mm to 420 mm, tolerance $\pm 1\%$
- Screw thread M85x1, except 4401-527-000-21 = M76x1
- Transmission $\geq 97\%$
- Transmission $\geq 75\%$ at VIS-range
- Laser-damage threshold coating up to 6 J/cm^2 at 1064 nm, 10 ns, 100 Hz
- Includes interchangeable protective glasses

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LINOS F-Theta-Ronar Lenses 1064/1030-1080 nm

- Fused-silica and optical-glass designs
- Telecentric versions available
- Focal lengths ranging from 70 mm to 420 mm, tolerance $\pm 1\%$
- Screw thread M85x1, except 4401-261-000-21 = M76x1
- Transmission $\geq 96\%$ with good performance in VIS-range
- Laser-damage threshold coating up to 40 J/cm² at 1064 nm, 9 ns and up to 1.5 J/cm² at 1030 nm, 285 fs
- Version 4401-xxx-000-26 with low absorption coating ≤ 20 ppm at 1064 nm
- Includes interchangeable protective glasses

LINOS F-Theta-Ronar 1064/1030-1080 nm, Quarzglas LINOS F-Theta-Ronar 1064/1030-1080 nm, Fused Silica

Nominal-brennweite (mm)	Max. Scanfeld (mm ²)	Scanwinkel $\pm \Theta_{x,y}$ max. (°)	Strahlenbündel-Durchmesser begrenzt bei 1/e ² \varnothing_{beam} (mm)	Bildpunkt-Durchmesser 1/e ² \varnothing_{spot} (µm)	Spiegelabstände (m1/m2) (mm)	Arbeitsabstand (WD) (mm)	Schutzglas	Artikel-Nr.
Nominal focal length (mm)	Maximum scan field (mm ²)	Scan angle $\pm \theta_{x,y}$ max. (°)	Beam diameter truncated at 1/e ² \varnothing_{beam} (mm)	Spot diameter at 1/e ² \varnothing_{spot} (µm)	Mirror distances m1/m2 (mm)	Working distance WD (mm)	Protective glass	Part No.
70 telecentric	30 x 30	± 12.4	14	10	17/12	87.5	PG22	4401-551-000-21
100 telecentric	43 x 43	± 12.2	14	15	17/28	137	PG14	4401-561-000-21
167 telecentric	84 x 84	± 14.8	20	17	26/28	215.4	PG16	4401-513-000-21
255	170 x 170	± 19.2	10	50	13/30	317.4	PG14	4401-499-000-26
340	205 x 205	± 17.7	14	51	17/29	441.6	PG16	4401-546-000-26
420	254 x 254	± 17.3	14	60	26/24	510.9	PG14	4401-508-000-26

LINOS F-Theta-Ronar 1064/1030-1080 nm, Optisches Glas LINOS F-Theta-Ronar 1064/1030-1080 nm, Optical glass

Nominal-brennweite (mm)	Max. Scanfeld (mm ²)	Scanwinkel $\pm \Theta_{x,y}$ max. (°)	Strahlenbündel-Durchmesser begrenzt bei 1/e ² \varnothing_{beam} (mm)	Bildpunkt-Durchmesser 1/e ² \varnothing_{spot} (µm)	Spiegelabstände (m1/m2) (mm)	Arbeitsabstand (WD) (mm)	Schutzglas	Artikel-Nr.
Nominal focal length (mm)	Maximum scan field (mm ²)	Scan angle $\pm \theta_{x,y}$ max. (°)	Beam diameter truncated at 1/e ² \varnothing_{beam} (mm)	Spot diameter at 1/e ² \varnothing_{spot} (µm)	Mirror distances m1/m2 (mm)	Working distance WD (mm)	Protective glass	Part No.
100	62 x 62	± 17.7	12	16	16/12	97.7	PG2	4401-302-000-21
100 telecentric	54 x 54	± 15.6	15	13	20/32	126	PG6	4401-464-000-21
160	99 x 99	± 17.7	12	26	16/12	176.2	PG2	4401-301-000-21
163	115 x 115	± 20.2	10	32	13/24	185.9	PG5	4401-261-000-21
254	157 x 157	± 17.7	20	25	26/23	296.2	PG6	4401-288-000-20
330	217 x 217	± 18.7	16	40	18/24	387.6	PG6	4401-360-000-21
420	291 x 291	± 19.8	15	55	30/16	494.2	PG6	4401-350-000-21

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LINOS F-Theta-Ronar Lenses 1550/1900-2000 nm

- Fused-silica and optical glass designs
- Telecentric versions available
- Focal lengths ranging from 100 mm to 437 mm, tolerance $\pm 1\%$
- Screw thread M85x1
- Transmission $\geq 95\%$ at 1550 nm
- Transmission $\geq 96\%$ at 1940 nm
- Laser-damage threshold coating up to 20 J/cm^2 at 1064 nm, 12 ns, 100 Hz
- Includes interchangeable protective glasses

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Protective Glasses

- Optimum protection for the optical system
- Coated on both sides
- High transmission for the corresponding wavelength or wavelength range
- High laser-damage threshold
- Short delivery time

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
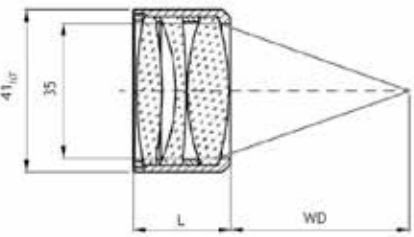
LINOS Focus-Ronar Lenses, 355 nm, Fused Silica

Focussing lenses are optimized for high precision applications, as used in laser systems for welding, cutting, drilling and structuring.

- Focal lengths ranging from 58 mm to 120 mm
- Flexible lens exchange without any adjustment
- Three-lens-element designs
- Full fused-silica designs
- Diffraction limited up to 17.5 mm ($1/e^2$) entrance beam diameter
- High clear aperture 35 mm
- Housing diameter 41mm
- Transmission $\geq 98\%$
- Damage threshold 5 J/cm² at 355 nm, 5 ns, 100 Hz

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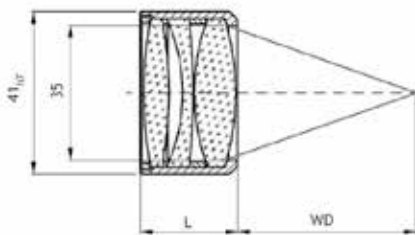LINOS Focus-Ronar Lenses 1064 nm/532 nm

Focussing lenses are optimized for high precision applications, as used in laser systems for welding, cutting, drilling and structuring.

- Focal lengths ranging from 58 mm to 120 mm
- Flexible lens exchange without any adjustment
- Three-lens-element designs
- High clear aperture 35 mm
- Housing diameter 41_{h7} mm
- The coating is optimized for 1064 nm, $T \geq 97\%$ and $T(532\text{ nm}) \geq 96\%$
- Good inspection performance at VIS wavelengths
- Laser-damage threshold 10 J/cm² at 1064 nm, 9 ns, 100 Hz
- Laser-damage threshold 6 J/cm² at 532 nm, 6 ns, 100 Hz

LINOS Fokus-Ronar 1064+532 nm, Optisches Glas LINOS Focus-Ronar 1064+532 nm, Optical glass

Brennweite (mm)	Länge L (mm)	Arbeitsabstand WD bei 1064 nm (mm)	Arbeitsabstand WD bei 532 nm (mm)	Part No.
Focal length (mm)	Length L (mm)	Working distance WD for 1064 nm (mm)	Working distance WD for 532 nm (mm)	Part No.
58	24.6	48.3	47.7	4401-505-000-20
77	18.9	72.2	71.5	4401-486-000-20
90	33.6	73.7	73.3	4401-490-000-20
120	24	110.7	109.8	4401-420-000-20

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Please look at www.qioptiq-shop.com for datasheets to our laser material lenses. The datasheets can be found at each product under Docs + Drawings. Our Laser Material Processing brochure is also ready to download here.

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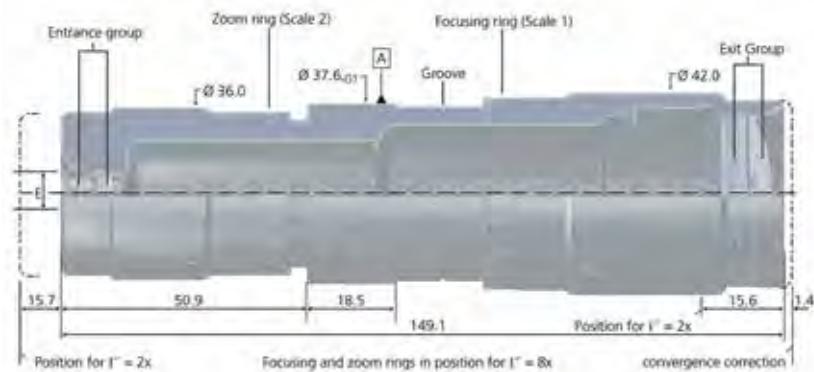
LINOS Variable Magnification Beam Expanders

- Continuous variation of magnification 2x...8x or 2x...10x
 - Choice between fused-silica or glass entrance lens
 - Continuous variation of exit-beam divergence
 - Wavelengths 355 nm, 532 nm, 633/780/830/980 nm or 1064 nm
 - Precise scales allow reliable settings and high repeatability
- Max. exit-beam diameter 31 mm
 - Max. length 165.5 mm
 - Max. diameter 42.0 mm
 - Mounting diameter 37.6_{-0.1} mm, reference on surface A

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LINOS Beam Expanders are optimally employed in conjunction with LINOS F-Theta-Ronar lenses for applications including:

- Laser structuring of foils
- Laser scribing of ceramic substrates
- Cutting of solar cells
- Micro drilling of sheet metal
- Marking of diverse materials with encodings

All LINOS beam expanders can also be implemented in reverse mode as beam reducers.

LINOS Variable Strahlaufweitungen 2x ... 8x
LINOS Variable Magnification Beam Expanders 2x...8x

Wellenlänge (nm)	Max. Eintrittsbündel-Ø bei 1/e ² Gauß-Strahl (mm) ¹⁾	Eintrittslinse aus	Part No.
<i>Wavelength (nm)</i>	<i>Max. entrance beam-Ø at 1/e² Gaussian beam (mm) ¹⁾</i>	<i>Entrance lens made of</i>	<i>Part No.</i>
355	3.4	Fused silica	4401-402-000-20
532	4	Fused silica	4401-446-000-20
532	8	Optical glass	4401-257-000-20
633/780/ 830/980	8	Optical glass	4401-258-000-20
1064	4	Fused silica	4401-359-000-20
1064	8	Optical glass	4401-256-000-20

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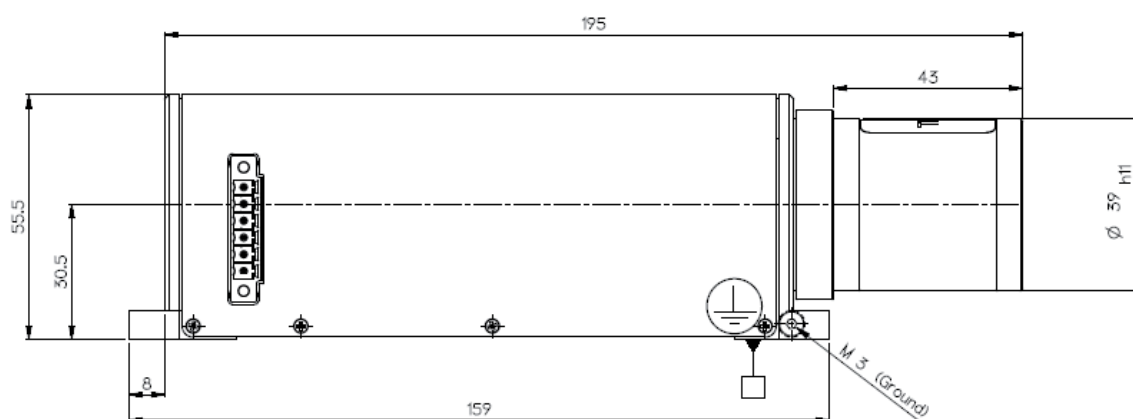
LINOS Motorized Beam Expanders

- Continuous variable magnification 2x...8x
- Fused-silica or optical-glass designs
- Continuous variation of exit-beam divergence
- Wavelengths 340-360 nm, 515-540 nm or 1030-1080 nm
- Software running on the Windows™ platform (XP, Win7, Win8)
- Reduce machine setup times by automatic change of magnification
- Maintains laser-protection class during readjustment of the beam expander
- All-in-one design, controller integrated
- CE and ROHS conform
- IP 20

- Exit-beam diameter: max. 31 mm
- 10 individual pre-sets for magnification and divergence
- Pointing stability < 0.5 mrad
- Fast adjustment from 2x to 8x about 7 s
- Mechanical dimensions: length 203 mm, width 58 mm, height 55.5 mm
- Mechanical interface via high-precision holes 6^{H7} (recommended) or mounting diameter 39_{H11}
- Different electronic interfaces: SubD9/RS232, USB 2.0, Phoenix Contact/RS232
- Baud rate: 9600 bit/s
- Power input 7–12 V, Phoenix Contact 7-24 V
- Entrance-beam diameter max. = 31 mm / zoom factor

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

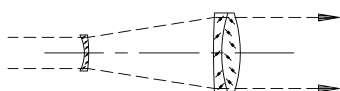
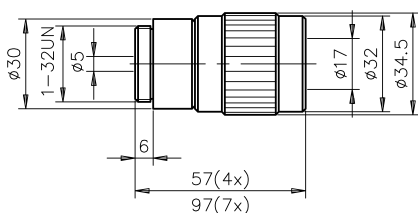
Laser Beam Expander System 4x and 7x

- Broadband anti-reflection coated with ARB2 for $\lambda = 450-700 \text{ nm}$
- To generate plane wave fronts
- To focus laser beams at long distances
- To reduce laser beam divergence
- To use in alignment work or optoelectronic control systems
- Equipped with focusable exit optics
- Fits directly to most lasers
- Adaptable to Microbench

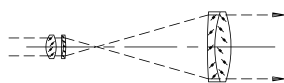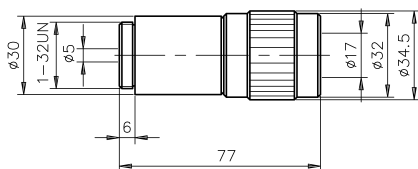- Entrance aperture: $\varnothing 4 \text{ mm}$
- Exit aperture: $\varnothing 16 \text{ mm}$
- Entrance element: plano-concave singlet
- Exit element: focusable from 0.5 m to ∞ (4x); focusable from 1.5 m to ∞ (7x)
- Mating thread: 1" x 1/32"

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Laser Beam Expander System 10x

- Broadband anti-reflection coated with ARB2 for $\lambda = 450-700 \text{ nm}$
- To generate plane wave fronts
- To focus laser beams at great distances
- To reduce laser beam divergence
- To use in alignment work or optoelectronic control systems
- Equipped with focusable exit optics
- Fits directly to most lasers
- Adaptable to Microbench

- Entrance aperture: $\varnothing 3 \text{ mm}$
- Exit aperture: $\varnothing 17 \text{ mm}$
- Entrance element: air-spaced achromat
- Exit element: focusable from 1.5 m to ∞
- Mating thread: 1" x 1/32"

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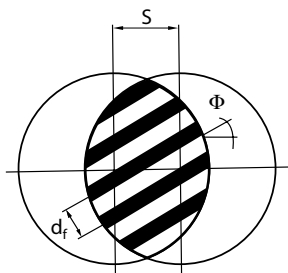
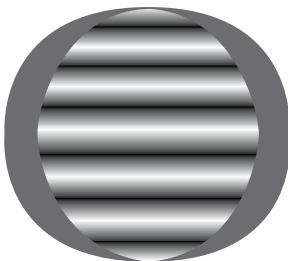
Please look at chapter Microbench, section Adapter regarding these accessories to our Laser beam Expander Systems.

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Shearcubes

The laser beam is being reflected at both front and back side of the wedge-shaped beam splitter plate. Both reflected beams are overlapped. The area of overlap is imaged on the dispersion plate of the Shear Plate Interferometer and can be observed easily. The laser beam is collimated, once the interference fringes are parallel to a reference line on the dispersion plate. If the beam expander is focussing or defocussing, respectively, the fringes are rotated clockwise or counter clockwise, respectively.

A quantitative analysis of the interferogram allows for the determination of the radius of curvature R of the wavefront. The laser divergence DIV can be determined easily, if radius of curvature R and diameter of the laser beam d_{Strahl} are known.

- For fast determination of laser collimation

Other applications:

- Determination of refractive index of plane plates (accuracy $\pm 1 \cdot 10^{-3}$)
- Determination of focal lengths of lenses (accuracy $\pm 3 \%$)
- Determination of radius of curvature of concave and convex mirrors (accuracy $\pm 3 \%$)
- Three versions available for beam diameters between 1 mm and 25 mm
- Compatible with LINOS Microbench, Nanobench, and Tube C systems
- Beam entrance height: 20 mm
- Robust and easy to use

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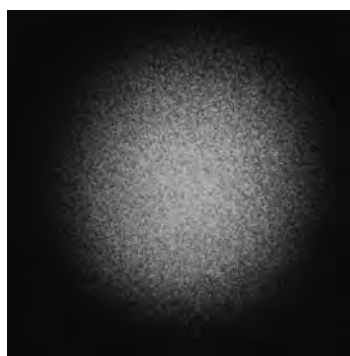
Laser Speckle Reducer


Speckle noise from a laser-based system is reduced by dynamically diffusing the laser beam. Electro-active polymers are used to oscillate a diffuser in its plane (x- and y-direction). This innovative principle is very power-efficient and completely free of mechanics offering years of lifetime. The LSR-3000 series integrates fully certified drive electronics powered through a single micro-USB connector. The LSR is available in two versions: the LSR-3005 and the LSR-3010 that exhibit a clear aperture of 5 mm and 10 mm, respectively.

You can find more information and datasheets at www.optotune.com


- Power supply: 5 V DC (micro-USB interface)
- Power consumption: 310 mW

CCD image of a laser spot without LSR-3005 on

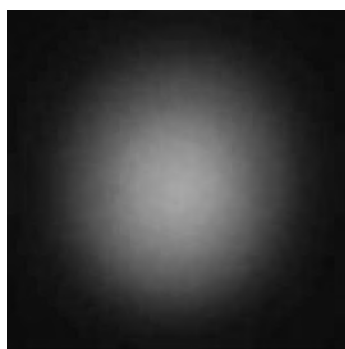CCD image of a laser spot with LSR-3005 on

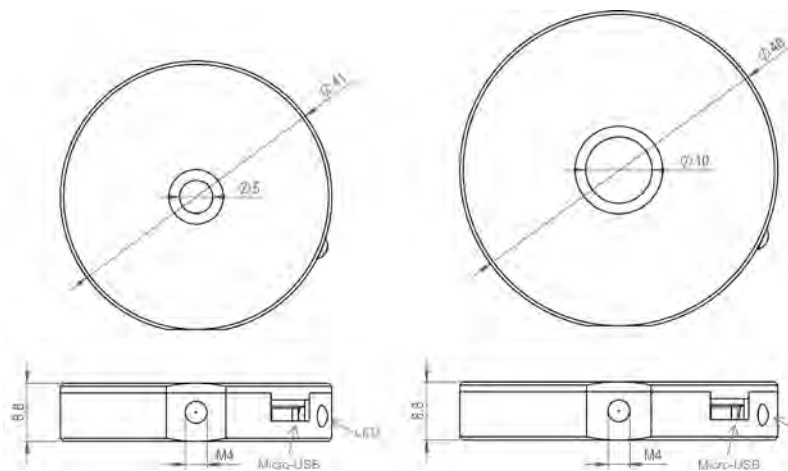Mechanical drawing of LSR-3005 (left) and LSR-3010 (right)

Laser Speckle Reducer LSR3000 Laser Speckle Reducer LSR3000

Produkt-bezeichnung	Durchmesser (mm)	Wellenlängenbereich (nm)	Freie Öffnung Ø (mm)	Streuwinkel (°)	Durchmesser (mm)	Dicke (mm)	Oszillationsfrequenz (Hz)	Part No.
Item Title	Diameter (mm)	Wavelength range (nm)	Clear aperture Ø (mm)	Diffusion angle (°)	Diameter (mm)	Thickness (mm)	Oscillation frequency (Hz)	Part No.
LSR-3005-6D-VIS	41	400-700	5	6	41	8.8	300	G399557000
LSR-3005-12D-VIS	41	400-700	5	12	41	8.8	300	G399558000
LSR-3005-24D-VIS	41	400-700	5	24	41	8.8	300	G399559000
LSR-3005-17S-VIS	41	400-700	5	17	41	8.8	300	G399560000
LSR-3005-6D-NIR	41	700 - 1100	5	6	41	8.8	300	G399561000
LSR-3005-12D-NIR	41	700 - 1100	5	12	41	8.8	300	G399562000
LSR-3010-6D-VIS	48	400-700	10	6	48	8.8	180	G399563000
LSR-3010-12D-VIS	48	400-700	10	12	48	8.8	180	G399564000

The LINOS Machine Vision Solutions

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Our quality criteria:

- *Excellent image quality over the entire image field*
- *Low distortion*
- *Rugged quality for use in industry*
- *High color correction in the visible range*

Machine Vision Solutions

Qioptiq manufactures lenses for all of today's image processing tasks. A broad range of high-quality standard lenses means LINOS lenses can be used in practically all conceivable applications in image processing.

With well-established expertise and many years of experience, our research and development department is ready to implement customer-specific production, from OEM solutions to finished modules, at any time. Contact us!

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Special Features:

Thanks to a sophisticated system of focusing aids and adapters, LINOS lenses can be connected to any camera on the market. Customer-specific interfaces can be integrated as well.

Ideal Areas:

Traffic supervision, wafer inspection, PCB and FPD inspection, packaging inspection, tablet control, foreign matter detection, bar code scanning, measurement of overhead lines.

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- Highest optical performance
- Large image circle up to 1 inch
- For pixel size even below 2 µm
- High numerical aperture

- Focal length: 12 ... 50 mm
- Magnification range: 0 ... -0.1
- Spectral range: 450-950 nm
- Focusing: manual, lockable
- Iris diaphragm: manual, lockable
- Filter thread: M35.5x0.5
- Lens diameter: 42 mm

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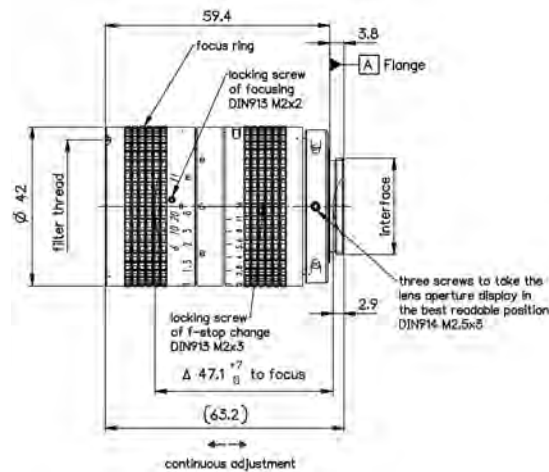
MeVis-C

Machine Vision lenses for Multi-Megapixel cameras

The LINOS MeVis-C lenses are specifically developed to be used with the highest resolution sensors available on the market. Exceptionally high resolution across the entire sensor ensures optimal performance for the most demanding applications.

High resolution is paired with low distortion, minimal light fall off and excellent chromatic correction across the entire spectral range of 450-950 nm. Premium glass and specialized coatings further combine to deliver very good color rendition. These lenses can resolve up to 200 lp/mm, even to the extreme corners of a 1" sensor. Only Mevis-C lenses combine all these positive optical characteristics and represent the perfect match for modern high-resolution sensors with up to 1" diameter.

The housing features locking screws for focus and aperture that prevent changes due to accidental contact or vibration and shock in harsh environments. Focus and aperture ring operate as smoothly as you would expect from a high-end lens and enable exact focusing and iris adjustment.

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MeVis-CF

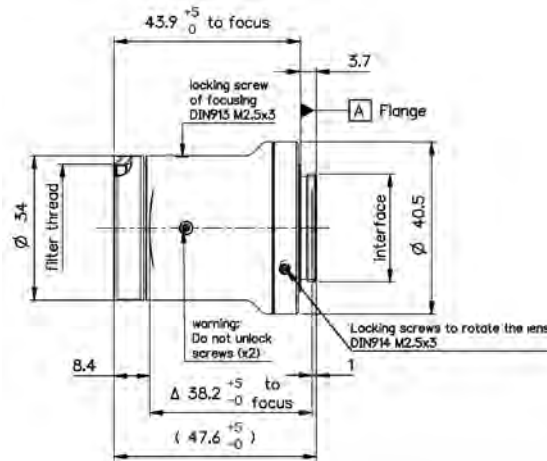
With the new MeVis-CF series, Qioptiq introduces a new line of lenses to serve applications requiring the best possible imaging quality from a compact, robust lens package. The optical performance is equal to the renowned MeVis-C series, but the mechanics are designed to be as simple as possible.

The major difference to the standard MeVis-C Series is the fixed aperture. This enhances durability, reduces size for compact installation and offers a more cost-effective design. The lenses are available in several versions offering a variety of apertures.

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- Highest optical performance
- Fixed aperture
- Large image circle up to 1 inch
- For pixel size even below 2 µm
- High numerical aperture
- Compact and robust mechanics

- Focal length: 16 ... 50 mm
- Magnification range: 0...-0.1
- Spectral range: 450-950 nm
- Focusing: manual, lockable

MeVis-CF MeVis-CF

Produktbezeichnung	Brennweite (mm)	Blendenzahl	Maßstabsbereich	Bildfelddurchmesser (mm)	Anschluss	Part No.
Item Title	Focal length (mm)	F-number	Magnification range	Image field of view (diameter) (mm)	Connector	Part No.
MeVis-CF 1.6/16	16	1.6	-0.1 ... 0	11	C-Mount	0020-009-000-20
MeVis-CF 2.8/16	16	2.8	-0.1 ... 0	11	C-Mount	0020-009-000-21
MeVis-CF 4.0/16	16	4.0	-0.1 ... 0	11	C-Mount	0020-009-000-22
MeVis-CF 25 mm f/1.6	25	1.6	-0.1 ... 0	16	C-Mount	0020-007-000-20
MeVis-CF 25 mm f/4	25	4	-0.1 ... 0	16	C-Mount	0020-007-000-21
MeVis-CF 25 mm f/2.8	25	2.8	-0.1 ... 0	16	C-Mount	0020-007-000-22
MeVis-C 35 mm f/1.6	35	1.6	-0.1 ... 0	16	C-Mount	0020-008-000-20
MeVis-C 35 mm f/2.8	35	2.8	-0.1 ... 0	16	C-Mount	0020-008-000-21
MeVis-C 35 mm f/4.0	35	4.0	-0.1 ... 0	16	C-Mount	0020-008-000-22
MeVis-CF 1.8/50	50	1.8	-0.075 ... 0	16	C-Mount	0020-010-000-20
MeVis-CF 2.8/50	50	2.8	-0.075 ... 0	16	C-Mount	0020-010-000-21
MeVis-CF 4.0/50	50	4.0	-0.075 ... 0	16	C-Mount	0020-010-000-22

NEW

MeVis-C/CF Traffic

High Resolution Inspection Lens with optimal VIS and NIR Performance

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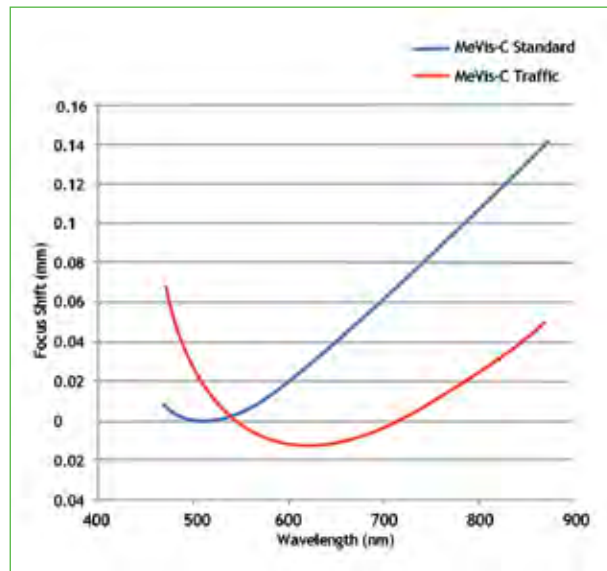
MeVis-C/CF Traffic

High Resolution Inspection Lens with optimal VIS and NIR Performance

The MeVis-C/CF traffic lenses feature optimized color correction to eliminate focus-shift between daylight and NIR illumination around 850 nm. This optimized optical design removes the need for refocusing when switching between wavelengths as the focal plane for both illumination cases is identical. The field-proven and well known mechanics of the MeVis-C and MeVis-CF lenses remain unchanged to ensure a robust lens ready for the most demanding industrial use cases.

- Optimized for visual spectrum and NIR with identical focus plane over broad wavelength range
- Highest optical performance
- Large image circle up to 1 inch
- For pixel size even below 2 μm
- High numerical aperture
- Spectral range: 450-950 nm
- Focusing: manual, lockable
- Iris diaphragm: manual, lockable
- Filter thread: M35.5x0.5
- Lens diameter: 42 mm

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Inspec.x M


Lenses for high resolution linescan and areascan applications

The inspec.x M Series closes the gap between the outstanding C-Mount MeVis-C lenses and the large format inspec.x L lenses.

Equipped with lockable manual focus and manual iris these F-Mount lenses are the ideal choice for sensors up to 35 mm format when a fast f/# is required.

The 1.4/50 mm lens is available for visible spectral range or near-infrared. The NIR version lens features a coating for the range from 900 to 1350 nm, making it an ideal choice for applications like electro- and photoluminescence.

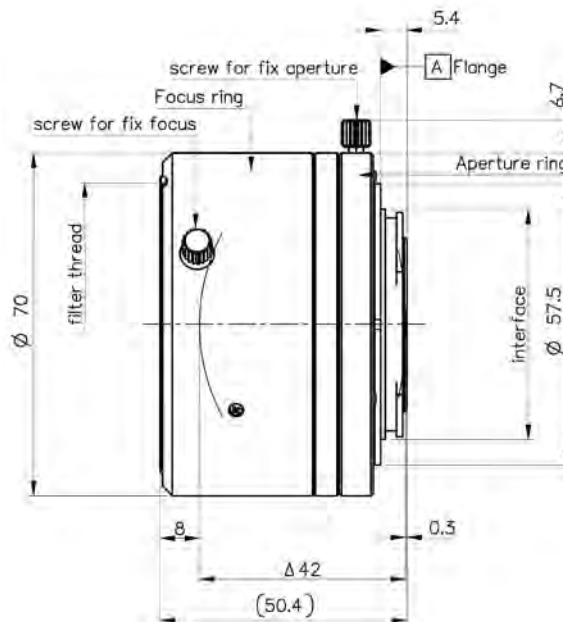- High resolution and contrast
- High magnification range
- High numerical aperture
- Magnification range: 0 ... -0.15
- Image circle: 43.3 mm
- Max. sensor size: 35 mm format
- Camera mount: F-Mount
- Focusing: manual, lockable
- Iris diaphragm: manual, lockable

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- Large image circle up to 82 mm
- Perfect match with 5 µm pixel size
- Full-metal barrel
- Lockable aperture

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Inspec.x L 5.6/105

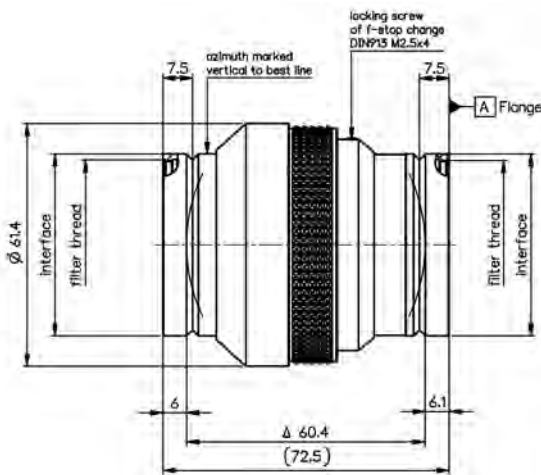
High-Resolution Lenses for Line-Scan Applications

The Inspec.x L Series was developed to meet the highest requirements in industrial image processing with very large sensors. This Lens Series shows even contrast and resolution over an image circle of up to 82 mm. The field-proven performance makes these lenses a perfect match for sensors like the popular 12k/5 µm and 16k/5 µm line-scan sensors. Very large area scan sensors also benefit from the high resolution, low distortion and excellent color correction of these lenses.

The four variants of the Inspec.x L 5.6/105 are optimized for magnifications of 0.33x, 0.5x, 0.76x and 1x respectively. The V-Groove interface at both ends of the lens enables adjustment to the best azimuth position and use of the lens in retro position for magnifications up to 3x.

All lenses feature lockable iris and are engineered for the use in harsh industrial environment.

- Focal length: 105 mm
- Aperture: 5.6 ... 11
- Working distance: 100-420 mm
- Spectral range: 400-750 nm
- Distortion: < 0.4%
- Camera mount: V-groove
- Compatible with all Inspec.x L series accessories

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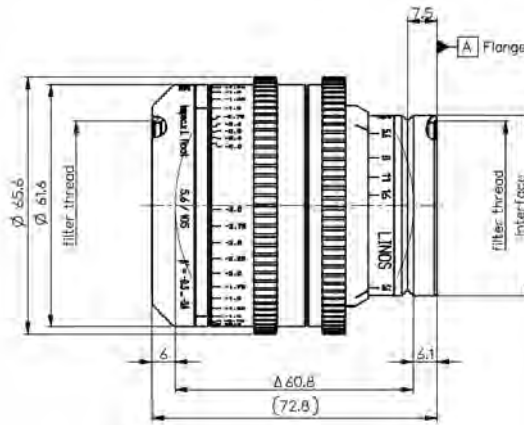
inspec.x L 5.6/105 float
High-Resolution Lens with Wide Magnification Range

The inspec.x L 5.6/105 float lens combines the field-proven imaging performance of the inspec.x L 5.6/105 Series with newly developed mechanics to cover a large magnification range with only one lens.

By using floating element optics the inspec.x L 5.6/105 float achieves high performance over an extremely wide magnification range from 0.3x to 3.0x. Full-metal housing makes the lens suitable for industrial use and geared rings facilitate the use in automated environments.

- Ideal for 12k and 16k sensors
- Geared rings for focus and aperture
- Large image circle up to 82 mm

- Focal length: 105 mm
- Aperture: 5.6 ... 11
- Extreme magnification range of -0.3x ... 3x
- Floating element optics
- Perfect match with 5 µm pixel size
- Working distance: 100-420 mm
- Spectral range: 400-750 nm
- Distortion: < 0.4%

inspec.x L 5.6/105 float
inspec.x L 5.6/105 float

Produktbezeichnung	Brennweite (mm)	Blendenzahl	Maßstabsbereich	Bildfelddurchmesser (mm)	Anschluss	Part No.
Item Title	Focal length (mm)	F-number	Magnification range	Image field of view (diameter) (mm)	Connector	Part No.
inspec.x L 5.6/105 float	105	5.6	-0.3 ... 3.0	82	V-groove	0703-114-000-20

- Herausragende MTF-Performance über den kompletten Bildkreis
- Beugungsbegrenztes Optikdesign
- Hohe numerische Apertur
- Outstanding MTF performance over the entire image circle
- Diffraction-limited optical design
- High numerical aperture

Inspec.x L 4/105

High-Resolution Lenses for High Magnification Imaging

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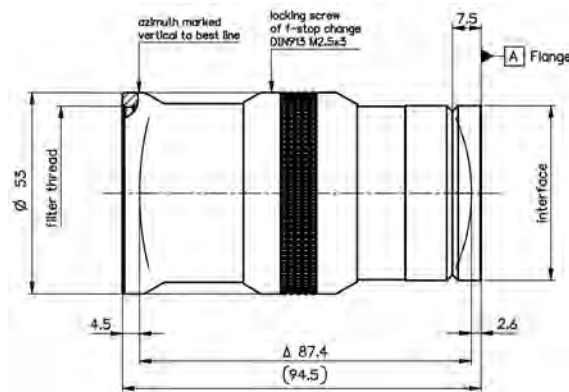
Inspec.x L 4/105

High-Resolution Lenses for High Magnification Imaging

The Inspec.x L 4.0/105 Series lens was developed to achieve an object resolution down to $2\ \mu\text{m}$. These lenses are designed for sensors with up to 82 mm length and achieve an enormous resolution of up to 200 lp/mm in the object plane. This high imaging quality is achieved across the complete image circle. Another advantage is the extremely small distortion, which is well under $1\ \mu\text{m}$ over the complete field of view.

The Inspec.x L 4.0/105 Series is the perfect match for modern, high-resolution $12\text{k}/5\ \mu\text{m}$ and $16\text{k}/5\ \mu\text{m}$ cameras for applications that require extremely high resolution. Three different magnifications of 3x, 3.5x and 5x are available. All three magnifications can also be ordered in a version this is optimized for a prism beamsplitter on the object side for coaxial illumination.

- Focal length: 105 mm
- Magnifications: 3x, 3.5x, 5x
- Image circle: up to 82 mm
- Spectral range: 400-750 nm
- Iris diaphragm: manual, lockable
- Mount: V-groove
- Wide range of mechanical accessories

- 3.7 ... -3.3
- 5.2 ... -4.8
- 3.3 ... -2.8
- 3.3 ... -2.8
- 4.0 ... -3.3
- 5.2 ... -4.8

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- Inspec.x L 4/105 -3x prism
- Inspec.x L 4/105 -3.5x prism
- Inspec.x L 3.5/105 -5x prism

Inspec.x L Prism Module

Beam Splitter Prism for High-Resolution Lenses

Coaxial illumination is very common for high-resolution applications. Qioptiq offers an easy-to-use prism module that is optimized for the use with the prism type Inspec.x L lenses. This prism module has a universal mount for fiber guides that are widely used for line-scan illumination. The setup of this system is very straight forward and results in a complete optical system of lens and illumination unit that achieves diffraction limited performance.

The prism module is designed for the lenses:

- Inspec.x L 4/105 -3x prism
- Inspec.x L 4/105 -3.5x prism
- Inspec.x L 3.5/105 -5x prism

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LINOS MachVis software is designed to help you find the right lens for your machine vision application.

- Starting with 4 simple parameters:
- Object size
 - Working distance
 - Sensor size
 - Camera mount

MachVis calculates the necessary optical parameters and selects all suitable machine vision components that meet the specifications.

For download and further information, please go to www.qioptiq.com/request-machvis as well as in chapter Optics Software.

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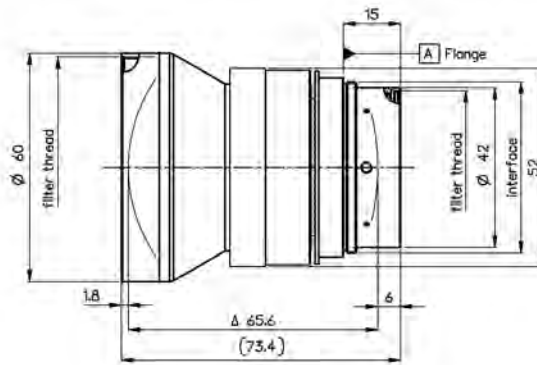
Inspec.x L 4/60 and 4/100

High-Resolution Lenses for Small Magnifications

The Inspec.x L 4/60 and 4/100 are optimized for the use from infinity down through magnifications of around 0.1x. In this range these lenses show exceptional contrast over a large sensor size of up to 70 mm. High contrast goes along with very good color correction and low distortion. Especially the 60 mm lens provides unusually high performance for such a short focal length and enables imaging of large objects in space constrained environments with large sensors.

Both lenses feature a threaded interface to use them with the popular Modular Focus helical mount that provides access to virtually all existing cameras via different camera adapters.

- High contrast for small magnifications
- Excellent color correction
- Focal length: 60 and 100 mm
- Magnification range: 0 ... 0.2
- Image circle: up to 70 mm
- Spectral range: 400-750 nm
- Iris diaphragm: manual, click-stop
- Mount: compatible to Modular Focus
- Wide range of mechanical accessories

Inspec.x L 4/60 und 4/100 Inspec.x L 4/60 and 4/100

Produktbezeichnung	Brennweite (mm)	Blendenzahl	Maßstabsbereich	Bildfelddurchmesser (mm)	Anschluss	Part No.
Item Title	Focal length (mm)	F-number	Magnification range	Image field of view (diameter) (mm)	Connector	Part No.
Inspec.x L 4.0/60	60	4.0	0 ... -0.2	70.8	M45x0.75	0019-002-000-49
Inspec.x L 4.0/100	100	4.0	0 ... -0.2	70.4	M45x0.75	0019-003-000-49

Singles
Achromats
Plane Optics
Mirrors
Thin Films
Coatings
Polarization Optics
Laser Optics

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Apo-Rodagon HR

High resolution lenses for line scan applications

The new Apo-Rodagon-HR is designed to close the gap between the high-resolution lenses of the inspec.x L Series, and the Apo-Rodagon-D lenses, both in terms of imaging quality and price.

The robust mechanics of this lens make it suitable for applications in the harshest environments. Thanks to the use of a fixed aperture there are no moving parts to shake out of adjustment due to vibrations in extreme applications. The standard aperture of 5.6 is the value at which the lens achieves the optimum performance. Other apertures are available on request.

The high resolution of 80 lp/mm at the sensor side in combination with the large image circle of 62 mm makes the Apo-Rodagon-HR 0.5x a very good match with the popular 12k / 5 µm line-scan cameras.

The Apo-Rodagon-HR 0.5x features the same focal length as the Apo-Rodagon-D 4.5/75 2x and can therefore replace this lens if higher resolution is needed.

- Ideal for 12 k / 5 µm sensors
- Compact and robust mechanics
- High performance / price value
- Large image circle up to 62 mm
- Magnification range: -0.65 ... -0.35
- Camera mount: V-Groove
- Iris diaphragm: fix
- Filter thread: M37x0.75
- Lens length: 47 mm
- Lens diameter: 46 mm

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Rodagon F

Precision optics for F-mount

The Rodagon-F Series was developed by Qioptiq to adapt precision industrial-grade optics to an F-Mount camera, with a great price/performance ratio. Now, for the first time, users can integrate world re-knowned Rodenstock image quality in 40-60 mm focal lengths directly onto F-Mount cameras. A revolutionary design eliminates all moving parts to offer exceptionally robust performance. The smooth focusing is locked with a massive retaining ring that is fixed with additional screws.

The Rodagon-F lenses are available in different versions with fixed apertures. The fixed aperture prevents accidental misadjustment of the iris or slowly shifting aperture values due to vibrations. The image circle of these lenses is 44 mm and is therefore large enough for sensors with 35 mm format and the popular 41 mm line sensors.

- Suitable for line-scan cameras and large imaging sensors
- Large image circle up to 46 mm
- High numerical aperture
- Focal length: 40 ... 60 mm
- Magnification range: 0 ... -0.5
- Spectral range: 400 ... 750 nm
- Iris diaphragm: fix
- Mount: F-Mount
- Filter thread: M40.5x0.5
- Wide range of mechanical accessories

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Rodagon M42

Rodagon M42 series with integrated focusing

The Rodagon lenses are successfully implemented in many different machine vision applications for many years already. Now, the Rodagon family grows further and Qioptiq sets a new standard with the Rodagon M42 series in terms of price-performance-ratio.

The new Rodagon M42 series is available with 40, 50 or 60 mm focal length. The lenses feature integrated manual focusing that can be adjusted precisely and fixed by a locking ring and additional clamping screws. Even strong vibrations will not cause defocusing of the setup. The flange focal distance of the M42 mount is specifically optimized for industrial line scan cameras like UNIIQA, Linea or Racer made by leading camera manufacturers to cover the complete magnification range of typical machine vision applications. The fixed aperture of the Rodagon M42 lenses further enhances the robustness as no accidental change of aperture setting is possible. The result is a simple and robust as well as cost-effective system comprising of camera and lens without any moving parts.

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APO-Rodagon D

High Performance Lenses for Large Imaging Sensors and Line-Scan applications

The high resolution of the Apo-Rodagon-D lenses makes them an optimum solution for cameras with a pixel size down to 7 μm . The high resolution is accompanied by ultra-low distortions and negligible color fringing. The lenses are optimized for a magnification range from 1:2 to 2:1.

- Specially designed for scanning applications and large imaging sensors
- Optimized for 1:2 to 2:1 imaging
- Large image circle up to 102 mm
- Focal length: 75, 120 mm
- Magnification range: -0.33 ... -3.0
- Spectral range: 400-750 nm
- Iris diaphragm: manual, click-stop
- Mount: M39x1/26" (Leica)
- Filter thread: M40.5x0.5
- Wide range of mechanical accessories

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Apo-Rodagon-N

Measuring Lenses for Large Imaging Sensors

The Linos measuring lenses developed by Qioptiq feature the highest resolution, excellent contrast, minimum distortion and color neutrality.

Apo-Rodagon-N

Lenses produce crisp, sharp images to the extreme edges of the object. Apo-Rodagon-N Apochromatically corrected lens designed to meet the highest requirements in an especially broad magnification range.

- Suitable for line-scan cameras and large imaging sensors
- Large image circle up to 100 mm
- High numerical aperture
- Adapter available for all common camera interfaces
- Focal length: 50 ... 105 mm
- Magnification range: -0.05 ... -0.5
- Spectral range: 400-750 nm
- Iris diaphragm: manual, click-stop
- Mount: M39x1/26" (Leica)
- Filter thread: M40.5x0.5
- Wide range of mechanical accessories

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Rodagon, Rodagon-WA

Measuring Lenses for Large Imaging Sensors

The Linos measuring lenses developed by Qioptiq feature the highest resolution, excellent contrast, minimum distortion and color neutrality. They sharply reproduce images all the way to the very edges of the object.

Rodagon

High-performance lens featuring a highly consistent imaging quality and broad magnification range.

Rodagon-WA

High-performance lens with extended-field viewing angle and a large resulting image circle.

- Suitable for line-scan cameras and large imaging sensors
- Large image circle up to 105 mm
- High numerical aperture
- Adapter available for all common camera interfaces
- Focal length: 28...135 mm
- Magnification range: -0.03 ... -0.5
- Spectral range: 400-750 nm
- Iris diaphragm: manual, click-stop
- Mount: M39x1/26" (Leica)
- Wide range of mechanical accessories

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Rogonar-S

Measuring Lenses for Large Imaging Sensors

The Linos measuring lenses developed by Qioptiq feature the highest resolution, excellent contrast, minimum distortion and color neutrality. They sharply reproduce images all the way to the very edges of the object.

Rogonar-S

Inexpensive lens with good imaging performance. Optimal price / performance ratio

- Suitable for line-scan cameras and large imaging sensors
- Large image circle up to 82 mm
- High numerical aperture
- Adapter available for all common camera interfaces
- Focal length: 50 ... 90 mm
- Magnification range: -0.08 ... -0.5
- Spectral range: 400-750 nm
- Iris diaphragm: manual, click-stop
- Mount: M39x1/26" (Leica)
- Filter thread: M40.5x0.5

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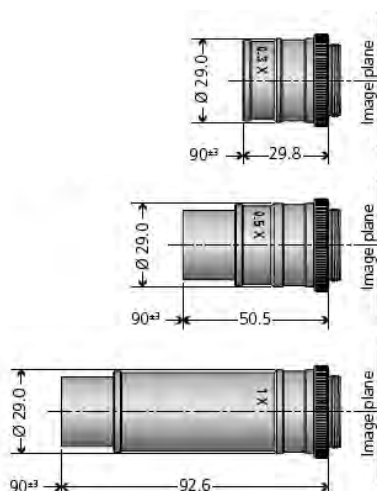
Macro CCD Lenses

Developed for C-Mount CCD Cameras

Macro Lenses developed by Rodenstock Präzisionsoptik for CCD Cameras, feature the highest resolution, excellent contrast, color neutrality and are virtually free of distortions. A large selection of Macro Lenses with magnification ratios from 1:3 to 4:1 is available which can be supplied with various fixed diaphragm aperture settings.

- High optical performance
- Huge range of magnifications available
- Designed for 1/2" and 2/3" sensors
- Image circle: 11 mm
- Lens diameter: 29 mm
- Working distance magnification 1:3 to 1:1: 90 mm
- Working distance magnification 2:1 to 4:1: 75 mm

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LINOS MachVis software is designed to help you find the right lens for your machine vision application.

Starting with 4 simple parameters:

- Object size • Working distance
- Sensor size • Camera mount

MachVis calculates the necessary optical parameters and selects all suitable machine vision components that meet the specifications.

For download and further information, please go to www.qioptiq.com/request-machvis as well as in chapter Optics Software.

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Compact Objectives

- Fixed aperture
- Compact design
- Optimized for an average image size
- Optimum image quality for object distance 700 mm
- Mount outer diameter: 16-25 mm

+ C-Mount Adapter für Kompaktobjektive finden Sie am Ende dieses Kapitels.
Please look at the end of this chapter regarding C-Mount adapter for compact objectives.

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FishEye Objective 190°



The lens was especially developed for use in pipe inspection. It delivers sharp images of tube walls, which allows software a two-dimensional "unfolding" of the tube. Damaged areas can now be more easily detected at a sewer inspection.

Due to the compact design, the objective can be universally adapted to cameras for the industrial sector. Therefore it is ideally suited for Machine Vision, Image Processing and Surveillance Systems.

- Developed for 1/3" sensor with total effective size 4.8 mm (H) x 3.6 mm (V)
- 190° diagonal FOV (incl. distortion)
- Outer diameter 20 mm
- Total length 19 mm
- Modifications on demand
- Focal length: 1.35 mm
- Diagonal FOV (incl. distortion): 190°
- Image circle: 3.3 mm
- F-number: f/2.8
- Object distance range through focusing: 60.0 mm - ∞
- Average transmission: ≥ 82 %
- Overall length (entrance surface to detector): 25 mm
- Back focal length (distance exit surface to detector): 5 mm

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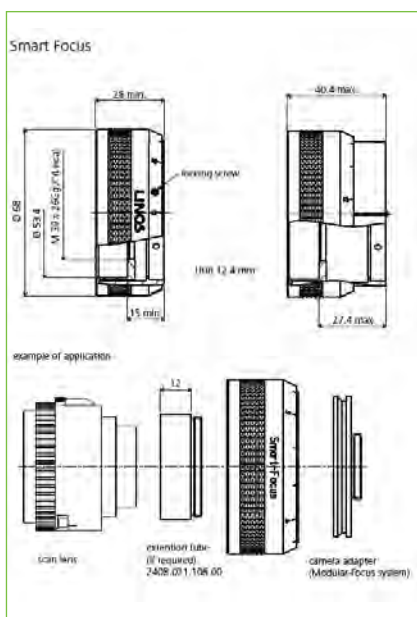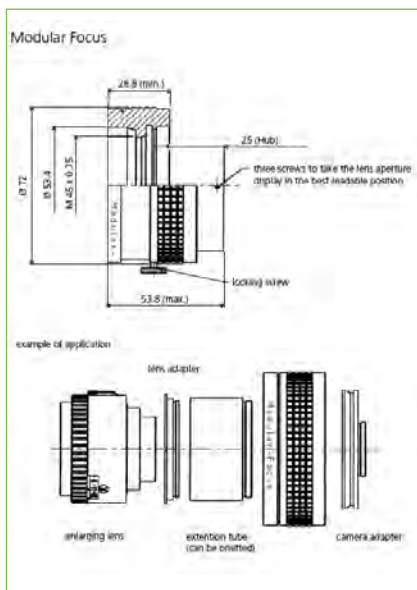Mechanical Accessories

Mechanical Adaptions

The LINOS Machine Vision Lenses have different threads and most have no focusing device. Therefore, we offer a complete range of mechanical accessories to adapt the lenses to all types of camera systems.

Modular Focus

A helical mount with a modular connection, can adapt the LINOS Lenses onto different camera threads. Using different lens and camera adapters, this modular application can connect various lens threads to various camera threads. The Modular Focus is designed to keep the lens from rotating during focusing.

Smart Focus

A low cost version of the Modular Focus with an M39x1/26" (Leica) lens thread. The diverse camera adapters of the modular focus are also used in this system. The difference between the Modular Focus and the Smart Focus is the shorter hub and that the lens is rotating while it is focused.

Focus Tube

The focus tube safely connects the inspec.x L lens-series to M72 and M95 camera mounts.

Modular Focus Helical mount

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Modular Focus Helical mount	2408-009-000-40
Camera adapter C-mount	2408-009-106-00
Camera adapter F-mount	2408-009-142-00
Camera adapter M42x1	2408-009-119-00
Camera adapter M48x0.75	2408-009-148-00
Camera adapter M58x0.75	2408-009-132-00
Camera adapter M72x0.75	2408-009-134-00
Camera adapter M90x1.0	2408-009-166-00
Camera adapter M95x1.0	2408-009-155-00
Extension tube 24.5 mm M45x0.75	2408-009-113-00
Extension tube 60 mm M45x0.75	2408-009-123-00
Extension tube 87.5 mm M45x0.75	2408-009-122-00
Extension tube 24 mm M72x0.75	2408-009-135-00
Extension tube 24 mm M90x1.0	2408-009-167-00
Extension tube 24 mm M95x1.0	2408-009-156-00
Lens adapter M32.5x0.5	2408-009-111-00
Lens adapter M39x1/26"	2408-009-118-00
Lens adapter M39x1/26"	2408-009-112-00
Lens adapter M45 V-Nut	2408-009-147-00
Retro ring M40.5x0.5 - M39x1/26"	2408-009-158-00
Retro ring M37x0.75 - M45x0.75	2408-009-152-00

Smart-Focus Helical Mount

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Smart focus	2408-021-000-43
Extension tube 12 mm M39x1/26"	2408-021-108-00
Extension tube 24 mm M39x1/26"	2408-021-109-00
Extension tube 48 mm M39x1/26"	2408-021-110-00
Extension tube 120 mm M39x1/26"	2408-021-111-00

Focus Tube inspec.x L 105

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
Focus tube M72 for inspec.x L 105 mm -0.33x and -0.5x	2408-012-000-31
Focus tube M72 for inspec.x L 105 mm -0.76x and -1.0x	2408-012-000-30
Focus tube M72 for inspec.x L 105 mm -3.0x and -3.5x	2408-012-000-47
Focus tube M72 for inspec.x L 105 mm -3.5x	2408-012-000-32
Focus tube M72 for inspec.x L 105 mm -5.0x	2408-012-000-33
Focus tube M95 for inspec.x L 105 mm -0.33x and -0.5x	2408-012-000-41
Focus tube M95 for inspec.x L 105 mm -0.76x and -1.0x	2408-012-000-43
Focus tube M95 for inspec.x L 105 mm -3.0x and -3.5x	2408-012-000-46
Focus tube M95 for inspec.x L 105 mm -5.0x	2408-012-000-45

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C-Mount Adapters for Compact Objectives

Mechanical Adaption

- Continuous adjustment via screws M2 (included in the scope of delivery)
- Thread length 4 mm
- Anodized

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Our quality criteria:

- *Precision engineering with innovative optical design*
- *Exact telecentricity on object and image side*
- *Diffraction limited optical performance over complete image circle*
- *Precise and dynamic AF*

LINOS Micro Imaging Solutions

For imaging of small objects most often traditional microscopes are being used. For integration into inspection systems and other machines these microscopes are often too complex and many features are integrated that are needed for machine integration. Qioptiq offers highly modular solutions for micro imaging that can be tailored to fit most applications without the bulk of a complete microscope.

This modularity is taken to the extreme with the Optem Fusion system that builds on the heritage of the Zoom 125 and Zoom 70XL families. With a multitude of configurations from fixed magnification to motorized 12.5:1 zoom including illumination and motorized focus this system is the right choice for demanding applications that require high imaging performance and ultimate flexibility.

With the mag.x system 125 Qioptiq offers an unparalleled optical system that is specifically designed to be used with modern imaging sensors of up to 57mm diagonal. mag.x system 125 offers high resolution already at low magnifications and supports object fields that are not possible with any other microscope optics.

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Special Features:

With the optional mag.x DIC module (differential interference contrast) even the most challenging inspection tasks can be solved.

Ideal Areas:

Inspection and processing of semiconductor, flat-panel display (TFT & OLED) and printed circuit boards (PCB), micro measurement & metrology, MEMS & nanotechnology, fluorescence microscopy, digital pathology and many more.

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mag.x system 125

High resolution imaging with wide field of view

The mag.x system 125 represents a new class of optical systems that enable microscope-like resolution with wide fields-of-view supporting modern high resolution sensors up to 57 mm diameter. It represents a completely new category of large field of view micro imaging lens systems. The variable system is fully modular and ideally suited for submicron-imaging in inspection and measurement applications.

The modular approach enables adaptation of the system to many applications ranging from a straight forward optical system without illumination to a full featured system including coaxial illumination and closed-loop autofocus.

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Tube Lenses

System magnification and maximum sensor size are the result of the combination of tube lens and objective lens. The current selection of exchangeable tube lenses (1x, 1.73x and 2.25x) allows the optimal use of line- and image-sensors with a diagonal of up to 25 mm, 43 mm and 57 mm.

All tube lenses are also telecentric on image side. Image space telecentricity helps to achieve exact measurement results and furthermore prevents brightness variations with sensors using micro-lenses.

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Base Units

Heart of the system is always the base unit. There are three different basic types of units available with additional options: base units with integrated illumination and optional turret, compact base units and base units with direct adapter.

All other components of the mag.x system 125 are attached to these base units, for example:

- Tube lens, objective lens and illumination tube
- Different polarizing filter sliders can be inserted into the base units
- Compact base units can be combined with a DIC-module in order to use the mag.x system 125 as a differential interference contrast microscope
- Base unit AF offers a universal port that is compatible to a variety of autofocus AF sensors, it can be combined with the optional piezo unit that enables through-the-lens AF

The base unit also provides the mounting interface of the whole system.

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LINOS MachVis software is designed to help you find the right lens for your machine vision application.

- Starting with 4 simple parameters:*
- Object size
 - Working distance
 - Sensor size
 - Camera mount

MachVis calculates the necessary optical parameters and selects all suitable machine vision components that meet the specifications.

For download and further information, please go to www.qioptiq.com/request-machvis as well as in chapter Optics Software.

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Piezo Units

High resolution systems, especially the used objective lenses, always feature small depth of field. Autofocus systems enable high resolution inspection of objects that are not perfectly flat.

The optional piezo unit is the active part for focusing. It can be combined either with the AF compact base unit that enables through-the-lens AF or with any other solution that detects the object position. The AF compact base unit offers a universal port that is compatible to a variety of AF sensors. This flexibility in autofocus and a perfectly homogeneous illumination allow operating a completely autonomous inspection system that can be controlled with a single control unit.

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Objective Lenses

The optical performance of the system is mainly defined by the objective lenses. These are the components that make the mag.x system 125 really unique. All of Qioptiq's highest end technologies are being used in manufacturing and testing these lenses.

mag.x system 125 stands out from conventional microscope systems by virtue of its low overall magnifications that transmit large object fields in high quality. This is achieved through numerical aperture (NA) values up to 0.32 far higher than those of other systems. High optical quality is not only ensured on the optical axis, but also is maintained over the entire sensor format.

Thanks to long working distances of up to a generous 24.8 mm and correction of the lenses without cover glasses, this system can be universally employed in technical inspection applications. The parfocal distance of the lenses is 125 mm and is origin of the name mag.x system 125.

Additionally exact object space telecentricity is precisely maintained, eliminating measurement errors due to 3-dimensional objects (height variations) or because of a variation in object position.

The optics is highly chromatically corrected in the spectral range of 430–700 nm. High contrast is maintained over the entire spectrum and no refocusing is required if the illumination wavelength is changed. Multispectral imaging becomes possible without any additional focus needs.

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Illumination

High-resolution inspection and measurement needs defined illumination that is optimally adjusted. This is why mag.x system 125 sports built-in coupled illumination for coaxial bright field illumination. The illumination system operates on the Koehler principle, with adjustable aperture and field stops. Thus the illumination and coherence parameters of the image (relationship of illumination and imaging aperture) can be appropriately adjusted for individual inspection tasks.

Additionally a special adapter is available for dark field illumination.

The coaxial bright field illumination is used in combination with the optional DIC-module in order to apply the mag.x system 125 as a differential interference contrast microscope.

Instead of applying the coaxial bright field illumination the system can be set up without integrated illumination as well using the direct adapter or customized solutions.

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Accessories

There are several camera adapters available. These mounts are used to connect the camera body of your choice to the tube lens of the mag.x system 125. The different female threads (M42, M72, M95) on the bottom are commonly used in the industry. The top side of the mount must match the camera body of the camera.

The interchangeable camera adapters allow the use of different format cameras, line scan, smart cameras, color or monochrome camera.

The mag.x system 125 can be combined with a DIC-module in order to use the system as a differential interference contrast microscope (to enhance the image contrast). It allows the inspection of transparent samples and objects with invisible features, where usual illumination techniques are not appropriate. For the DIC-operation one needs a compact base unit with two polarizing filters completed by an optional DIC modul. The handling and adjustment of the DIC-system is easy and user-friendly.

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Optem® FUSION System

Extreme Micro-Imaging Versatility

Optem® FUSION is the next generation of precision micro-imaging lens system. Modular design and infinity corrected optics combine for inspection, machine vision, monitoring and non-contact metrology applications. Designed for ultimate configuration flexibility to meet your exacting needs.

Features:

- Modular, micro-imaging lens system
- Change-out modules to modify form, function and performance
- User-configured for versatile 7:1 and 12.5:1 zoom imaging or economical fixed magnification – all in one system
- Optimized for most any camera format and mount type, including C, CS, F, 4/3 and other Bayonet mounts
- Programmable, touch-pad controller or board level electronics available to drive motorized options
- Improved optical design extends useable fields-of-view over previous Optem lens families

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Optem Fusion Basic Modules

The modular, micro-imaging lens system consists of the following types of components

- Camera adapter
- Camera tubes
- Zoom modules
- Lower function modules
- Lower lens and objectives

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Camera Mounts

Various Camera Mounts allow nearly all kinds of cameras to be mounted to the Camera Tubes that simply thread on to the T2 thread provided by the selected Camera Tube. On the other side standard connectors are available e.g. C/CS or F-mount to connect different cameras to the system.

- Compatible camera adapter allow the connection of different types of cameras

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Camera Tubes

FUSION Camera tubes provide the final magnification and image size for the optical system. They are adjustable to provide both par-focality and par-centricity to the image for best performance.

They are centerable with set screws or thumb-screws. The selection of tubes allows the best match of the camera to meet the application. Mini tubes make the whole system much more compact, in addition they show a reduces distortion.

- Common diameter through entire product line allows consistent mounting 1.625" (41.28 mm)
- Improved adjustment for parfocality and par-centricity
- Increased image sizes up to 43 mm diameter
- Large selection of different camera tubes
- Quick attachment and orientation
- Same tube used with right angle configurations
- Robust design
- Custom designed lenses for utmost optical quality and performance
- Perfectly matched for all FUSION combinations
- Same camera tubes are used for fixed and zoom configurations
- Preserves the numerical aperture of the lower modules
- Provide larger FOVs due to use of larger format cameras
- Supplemental magnifications from 0.4x to 3x
- Focal lengths from 80 mm to 600 mm
- Mini tubes designed for minimum length
- Transmission over the entire visible range

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Zoom Modules

There are two zoom ranges available with FUSION zoom core modules, both designed for the highest optical resolution and performance available. The 7:1 and the 12.5:1 zoom ranges were specifically designed for micro-inspection and imaging applications requiring both precision zoom quality and optical flexibility. Both zooms are available in three configurations; manual, detent and stepper motorized.

The zooms are uniquely designed for infinity on both ends for utmost configurability and ease of customization. When the zooms are used in any system, the magnification range is easily determined by the product of the lower lens magnification, the camera tube magnification and the reading on the zoom dial.

- Robust design
- Ease of attachment into system
- Customer convertible between 3 configurations (manual, detent, stepper)
- Conversion kits available
- Common diameter
- Meticulously designed for zoom repeatability, optical performance and constant focus throughout the zoom range
- Enclosed zoom gears

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Lower Function Modules

The vast selection of FUSION Lower Function Modules allows for the best solution to your needs. These modules are designed to allow integration of different lighting techniques, focus movement and lower lens options.

- Basic lowest cost mounting of the lower (taking) lens
- Internal focus modules
- Coax illumination and coax with focus modules
- Motorized versions
- All modules accept every lower lens or objective as well as attachment of ring lights
- Either cube accessory module attachment
- Fluorescent cube holder for large field applications

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Lower Lens

The FUSION lower lenses and objectives were all designed for seamless integration into the infinity optical portion of either zoom or fixed optical systems. There are considerable lower lens modules and microscope objectives to utilize in the FUSION system. This FUSION system does not "stack" additional auxiliary lenses into the system to vary the magnification and working distance, saving space and cost.

- Lower lens working distances from 32 mm to 490 mm
- Lower lens numerical apertures from 0.19 to 0.0036
- Objectives from 2x to 50x (require adapter)
- Objective numerical apertures from 0.055 to 0.6
- Easily changeable at any time
- Custom designed for best optical performance

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LWD Objectives

These objectives when attached to the FUSION system provide the utmost in resolution and working distance for demanding applications requiring higher magnification levels. The five M plan infinity corrected objectives provide flat field imaging for both zoom and fixed magnification needs.

These objectives are suitable for brightfield illumination as well as the specifically designed fiber optic ringlight for three dimensional objects. There are three HI resolution objectives for maximum resolution within the center portion of the image.

- Sufficient working distance from 13 to 34 mm
- Infinity corrected
- NA from 0.055 to 0.6
- Standard thread adapter required for mounting

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These LWD objectives are not just suitable for use with the Optem system. As they feature M26x36T threads, the objectives are also ideal for integration versatility with many metallurgical microscopes and industrial inspection systems wherever high resolution and long working distances are indispensable.

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Fetura+

The next generation Advanced Zoom Lens FETURA+ is a self-contained, programmable optical zoom lens that represents the apex of variable magnification performance and technology.

The most significant improvements over the first generation FETURA lens occur in the optical design. FETURA+ is infinity corrected on both sides of the zoom to permit compatibility with Qioptiq's Optem FUSION lens system accessories.

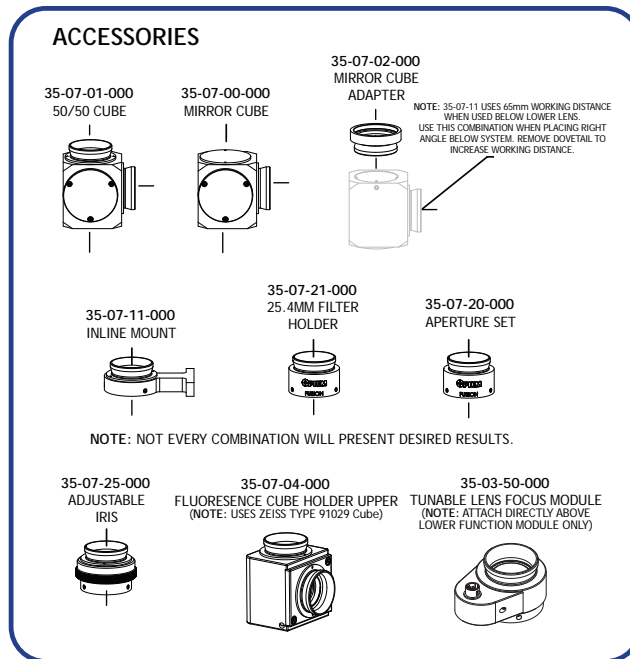
The FETURA+ optical redesign features improvements in optical performance and enhanced contrast, especially in the edges of the FoV.

Further improvements have been implemented in mechanics and electronics to achieve enhanced focus repeatability and mechanical stability for the most demanding applications and imaging environments. On the firmware side the continuous zoom feature was added, it enables zooming with a constantly focused object for a live view of the zooming progress – zoom speed can be adjusted to meet your viewing requirements. This can also be realized with a simple external potentiometer as well to achieve headless operation without a PC.

The unique linear-rail design and robust engineering ensure consistent repeatability and precision throughout an impressive 1-million zoom cycles.

- *Enhanced optical performance*
- *Continuous 12.5:1 optical zoom*
- *Improved repeatability*
- *Compatibility with all Optem FUSION accessories and components*
- *Adjustable optical zoom speed*
- *Extended service life - 1-mil cycles*

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Fokusvariables Linienmodul Tunable Lens Focus Module

Tunable Lens Focus Module

The addition of the shape changing polymer tunable lens Optotune EL16-40 to the FUSION product line provides many advantages for remote focusing.

The tunable lens module with a 16 mm aperture maintains good resolution with any of the FUSION systems from fixed magnification to either of the 7:1 or 12.5:1 zoom systems. The tunable lens is inserted in collimated space just above the lower lens or objective and changes the nominal working distance of a lower lens.

- Large focus range around the zero position
- Integrated temperature sensor to maintain good repeatability and focus on in-situ applications
- Operating temperature ranges from -20 to +65 °C
- Operating life exceeds 1 billion cycles
- Specially selected tunable lenses with reduced wavefront error compared to the standard version
- Including USB controller stick with software for easy installment
- Adds about 27 mm to the overall length of any system and eliminates the need for any lower focus module
- Spectral range of 450 nm to 950 nm provides an average of 95% transmission.
- Due to the large aperture, using at any angle other than vertical will produce strong coma in the image.

Produktbezeichnung	Part No.
Item Title	Part No.
Tunable lens focus module	35-03-50-000

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Cubes


Two versions of right angle cubes are available for the FUSION system, a mirror cube module and a 50/50 beamsplitter module. Both modules can be used for multi-imaging functions or space saving requirements.

- Both cubes can be attached to several locations
- 50/50 cube allows splitting of the system for many applications (dual Magnification and cameras eyepiece addition, etc.)
- Combine them for parallel dual cameras both with correct image orientation

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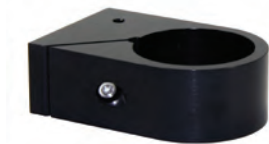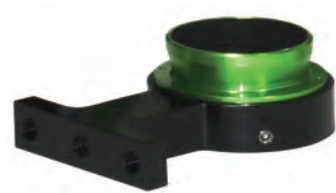
Mounts


The best mounting of the FUSION system requires rigid mounting of the optics. Several mounting accessories are available

- Camera tube mount anywhere along the tube
- Inline mount for between modules
- Both above provide a coplanar mounting surface with mounting holes
- Round "donut mounts" for installing in aftermarket microscope stands and stereo pods

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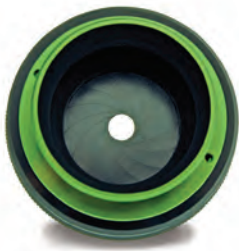Aperture Set, Filter Mount and Adjustable Iris Diaphragm

These modules provide customer adaptation at any time, the means to adjust the resolution, depth of field, intensity to camera or filtering. There are five apertures provided in the set and any 25 mm or 1 inch filter can be used.

- Used for adjusting intensity to the camera as well as numerical aperture and optical performance
- Filtering wavelengths to the camera
- Insert almost anywhere in the system

Einstellbare Filterbefestigung, and Apertur-Set
Adjustable Iris, Filter Holder and Aperture Set

Produktbezeichnung	Part No.
Item Title	Part No.
Adjustable iris	35-07-25-000
Filter holder 25.4 mm	35-07-21-000
Aperture set (required for fixed magnification systems)	35-07-20-000

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Fluorescent Cube Holder

This lower function module provides the means to use a standard Zeiss 91029 cube for large field fluorescent imaging and can be used with nearly any lower lens or objective.

- Easy integration of filters and dichroic mirror
- Less expense than a standard fluorescence microscope
- Easily adaptable for liquid fiber light guides for illumination excitation

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Infinity Macro Lens or Wide Field Lens


This attachment when combined with the 2x lower lens and a 15 mm focus module allows the customer to focus on objects from 50 mm to infinity. It provides a larger field of view (FOV) than the other lower lenses at a similar working distance.

- Easy to install
- Works with any fixed magnification or zoom configuration
- Quickly threads on the lower module

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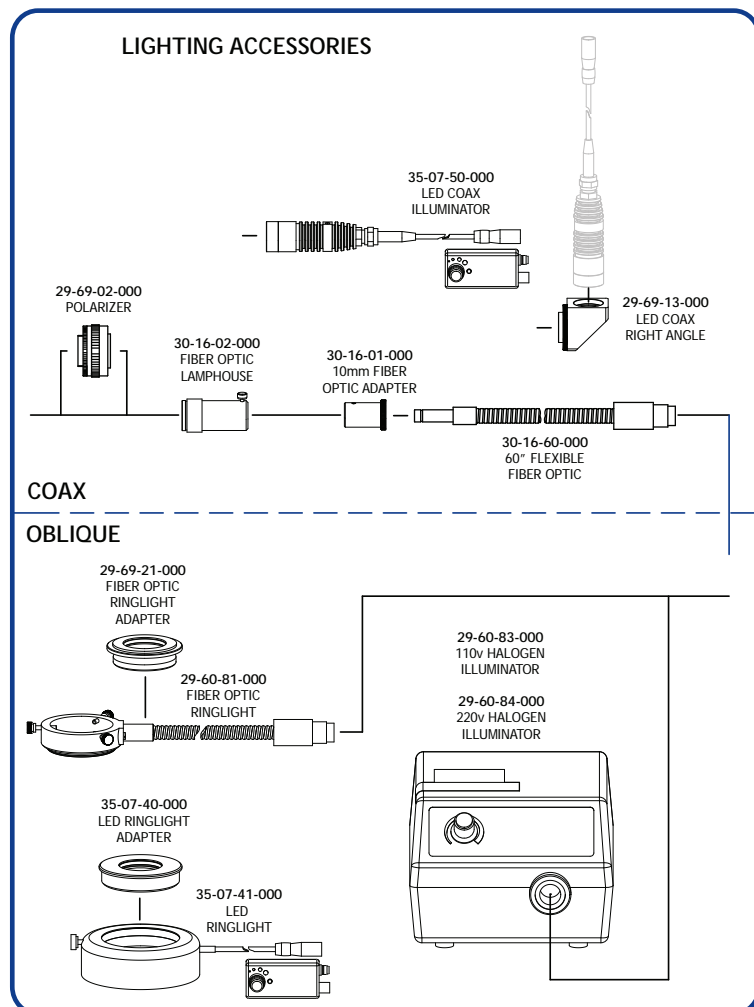Optem Fusion Beleuchtung Optem Fusion Illumination

Singlets
Achromats
Plane Optics
Mirrors
Thin Films
Coatings
Polarization Optics
Laser Optics
Machine Vision Solutions

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Optem Fusion Components for Illumination

FUSION illumination has been designed for easy integration into the product line. Several versions of both oblique (ringlight) and coaxial (brightfield) illumination are available. These are versatile for all kinds of imaging be it semiconductor, metrology, biotech, laboratory etc.. Our 150 watt halogen illuminator fits the flexible fiber as well as both fiber optic ringlights.

- Easy integration and mounting of components
- Minimal footprint
- Use of cool LED or fiber optic illumination
- Easily adaptable to third party illumination sources

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Optem Fusion Motorization

The newly designed stepper controller features programmable touchpad controls as well as a small clean footprint for space saving needs. For OEM users there is an OEM controller circuit board available for internal mounting to customer's equipment.

- These Controllers allow the control of two FUSION stepper motors be it zoom or focus or any combination thereof
- Stepping speed is user controllable
- 5 presets to automatically set predetermined zoom and focus settings
- 24 V DC required for OEM version

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A-Zoom2 10X Single-Objective Microscope

The A-Zoom2 10x single-objective zoom optical brightfield microscope delivers continuous 10:1 variable power magnification for microelectronics inspection, R&D imaging, testing and analysis.

Originally designed for electronic probing applications, A-Zoom2 features superior optics to locate and inspect the finest electronic traces using even finer electronic probes, A-Zoom2 features superior optics capable of resolving feature geometry in the sub-micron range.

- Manual 10:1 Zoom Optics
- Ergonomic Single-Objective design
- Select from a variety of Optem LWD Objectives
- Long-working distance maximize working space around subject
- Onboard CCD for video imaging
- 23 mm eyepieces with interocular adjustment
- Halogen coaxial illuminator with controller
- 50 mm heavy-duty focus block
- Easy integration with most prober stations
- Compatible with LINOS X95 Rail System and Stages for benchtop mounting
- Automated models available
- Laser-integrated models available

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Available for all A-Zoom Systems on request:

- Stable and rugged stand, as shown in the picture
- Motorized version
- Polarizer/Analyser combination
- More accessories

Please ask our specialist. Just send an email to sales.qioptiq@excelitas.com and he will contact you.

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A-Zoom2 40X Single-Objective Microscope

The A-Zoom2 40x single-objective zoom optical brightfield microscope delivers continuous 40:1 variable power magnification for microelectronics inspection, R&D imaging, testing and analysis.

Originally designed for electronic probing applications, A-Zoom2 features superior optics to locate and inspect the finest electronic traces using even finer electronic probes, A-Zoom2 features superior optics capable of resolving feature geometry in the sub-micron range.

- Manual 40:1 Zoom Optics
- Ergonomic Single-Objective design
- Select from a variety of Optem LWD Objectives
- Long-working distance maximize working space around subject
- Onboard CCD for video imaging
- 23 mm eyepieces with interocular adjustment
- Halogen coaxial illuminator with controller
- 50 mm heavy-duty focus block
- Easy integration with most prober stations
- Compatible with LINOS X95 Rail System and Stages for benchtop mounting.
- Automated models available
- Laser-Integrated models available

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A-Zoom Micro

The New A-Zoom Micro represents exceptional imaging value and economy for manual probe microscopy. A-Zoom Micro features a space efficient single-objective design which maximizes throughput, eliminates cumbersome nose turret manipulation and reduces risk of damage to fragile and high-value subjects.

Ideal for circuit board, semiconductor and flat-panel display probing, A-Zoom Micro is designed for streamlined integration onto most any probing station. It is also fully compatible for mounting to LINOS X95 Rail Systems and stages to provide a versatile high-magnification inspection system without sacrificing features, function or performance.

- Compact design
- Manual 7:1 parfocal optical zoom
- 2x Quick-slider extends magnification
- Ergonomic design
- 10x eyepieces
- Optimized for 1/2" C-Mount Cameras
- Easily configured for matching fields-of-view through eyepieces and monitor
- On-Board 10 W LED Coaxial illumination
- Heavy-duty focus block offers 50 mm fine-focus
- Optimized for use with a range of Optem LWD M-Plan APO Objectives

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Our quality criteria:

- *Extensive experience*
- *Top production facilities*
- *All sites benefit from ISO certifications*
- *Top production facilities*
- *Custom design support*

Our Capabilities

Qioptiq has top-qualified people and the state-of-the-art equipment to get the job done right. Qioptiq has project managers, R&D teams, sales advisors, and manufacturing and assembly capabilities in the United States, western Europe, eastern Europe and Asia. That means we can always find the right way, the right speed and the right price to respond to your needs.

We have over 100 years' experience in optics and photonics to offer our customers. Our close collaboration with universities and research facilities throughout the world, our efficient production processes and our worldwide customer base in multiple market sectors have enabled us to successfully set standards in the industry. Our renowned high quality is achieved thanks to our state-of-the-art production facilities and highly trained staff. Our expertise ranges from fast and cost-effective prototyping to consistent and reliable mass manufacturing — if necessary up to thousands of components over short periods of time.

With our highly skilled technical experts located on three continents, Qioptiq has all the skills required to provide you with fully customized system and design support. We are happy to apply our experience and our expertise to make your project successful.

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Special Features:

Two of our sites in Germany have extensive clean room technology

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Our Site in Asslar

<i>Micro Optics, General</i>	480
<i>Customized Micro Optical Components</i>	481
<i>Customized Opto-Mechanical Assemblies</i>	481
<i>Customized Micro Objectives</i>	481

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Our Site in Regen

Glass production has been at home in the Bavarian Forest for decades. This was just one more reason for Qioptiq to replace the aging plant with a new, state-of-the-art production site in 2008.

This plant is setting new standards with its versatility in production and the flexible deployment of their highly qualified skilled workers and engineers. The employees here on our two assembly lines are trained in a range of disciplines, and can be deployed wherever they are most needed.

At our site in Regen, 250 highly qualified staff are engaged in production using state-of-the-art manufacturing equipment

Starting with an individual work plan for each optical component, the combination of different manufacturing steps such as grinding, polishing, centring, quenching and tempering, precision bonding and edge varnishing create precision-optical elements to meet the highest requirements.

A glass technology laboratory for the determination of specific glass properties of the different materials is here to complement the production process.

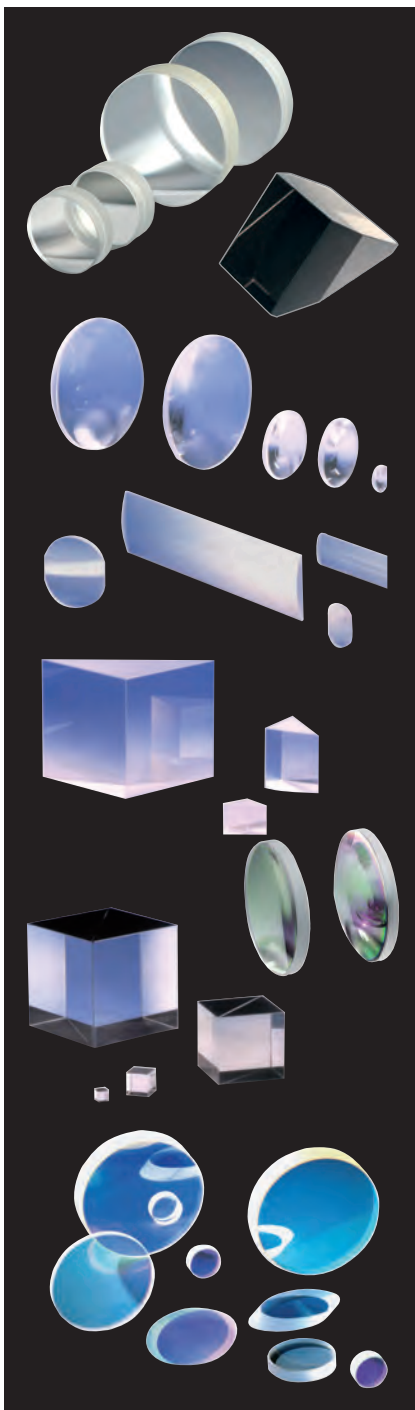
During assembly, complex optical systems for a wide range of applications are assembled from these materials together with electronic and mechanical components.

Precision Optical Components (POC) from Regen


Special production and series batches outside the catalogue range can be realised in our lens production at the Regen site.

We only require drawings with the geometries and optical specifications.

Technical advice on the applications is of course also available from one of our precision opticians. You need a special coating for example? Our coating designers will be happy to develop the optimized coating with you.

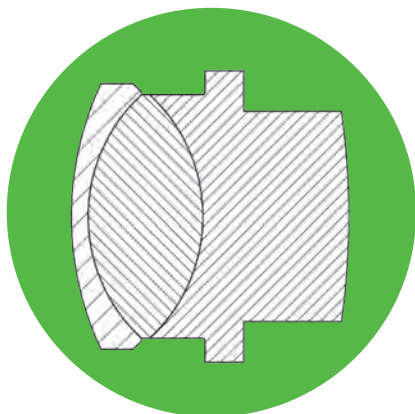
Singlets
Achromats
Plane Optics
Mirrors
Thin Films
Coatings
Polarization Optics
Laser Optics
Machine Vision
Solutions
Micro Imaging
Solutions

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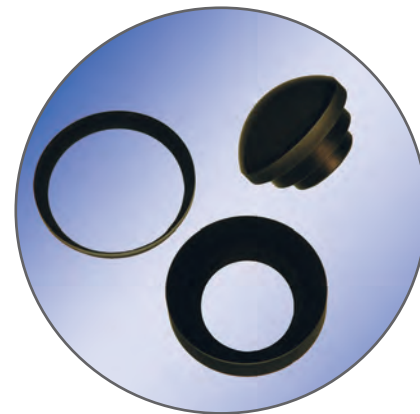Randlackieren / Edge varnishing

Capabilities

- Diameter: 3 to 400 mm
- Coating for all applications, new coating design if needed
- Special cooling for all glass materials
- Optical fine cement
- Documentation of all measured values
- Edge varnishing

Wolfgang Wilke, Factory Manager of our Regen Site

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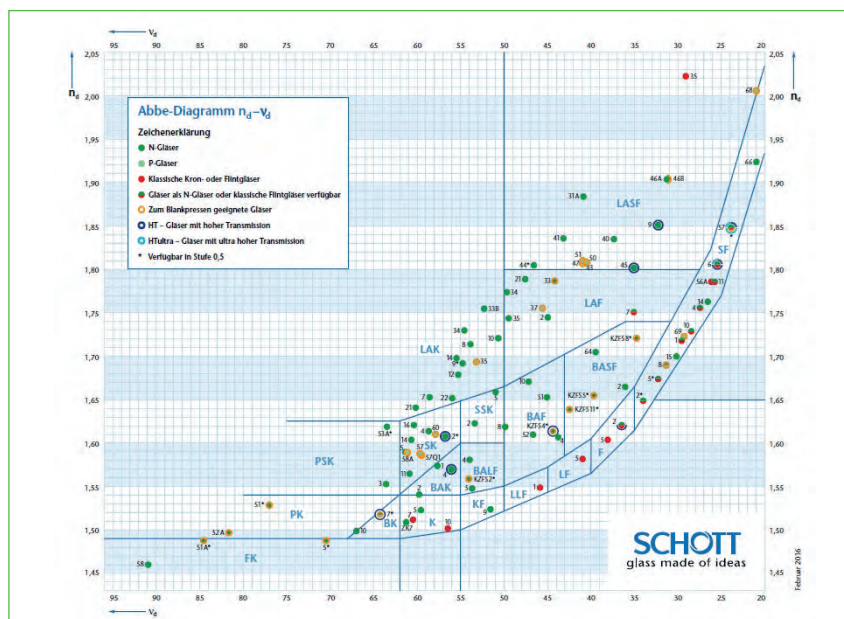
"Our mission is to fulfil the increasingly stringent requirements of geometrics, surface quality and cosmetics by (further) developing deterministic processes. Magnetorheological finishing and ion beam polishing is gradually becoming much more established. Welcome to the world of nanometers! Whereas in the past this degree of geometric precision was achieved almost entirely through the specific knowledge of polishers, today we are deploying ever more precision CNC machines and techniques. With the mastery of these newer processes, non-spherical geometrics – such as rotationally symmetric aspheres and freeform surfaces – are increasingly moving into the spotlight, including from an economical standpoint".

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Materials


- Complete portfolio of all well-established glass materials from SCHOTT and other premium glass suppliers like O'Hara
- Infrared materials (non-toxic), e.g. Germanium, Silicon
- Fused silica
- Sapphire
- Crystals (e.g. CaF_2 , MgF_2)

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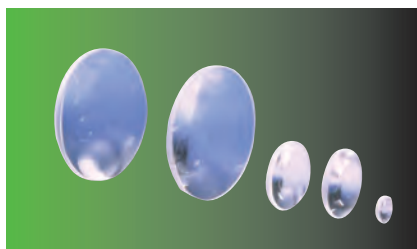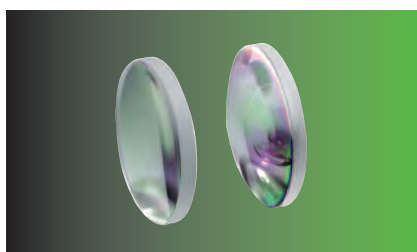
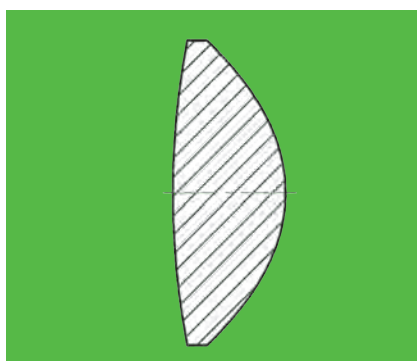
Peak-to-Valley Toleranzen Peak-to-Valley Tolerances

Diameter Optics up to (mm)	10	16	25	40	60	80
Peak-to-valley (µm)	5	5	15	30	70	70

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Optics

The following products can be produced for you:

Individual lenses (spheres)

- Diameter: Ø 3 to Ø 400 mm
- Form error: 3/0.1 (0.03)
Rmsl 0.002
- Cleanliness: to 5 / 0.025
- Diameter tolerance: to h6
- Thickness tolerance: to ±0.01

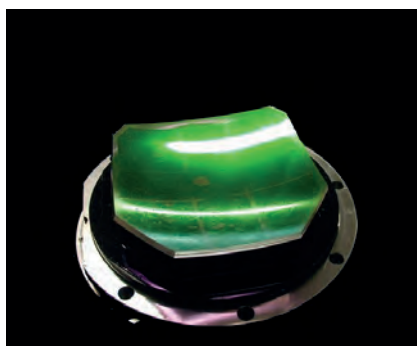
Aspheres, manufactured by conventional methods

- Grinded and polished
- Tolerances depending on the geometry

Aspheres, precision molded

- Glass types: B 270 and F2
- Possible diameter: Ø 3 mm up to Ø 80 mm

The difference between the maximum and minimum distance between two points is called peak-to-valley. In the case of optical glass, it denotes the mismatch between the actual or "test" surface and the ideal surface.

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
Freeform surfaces

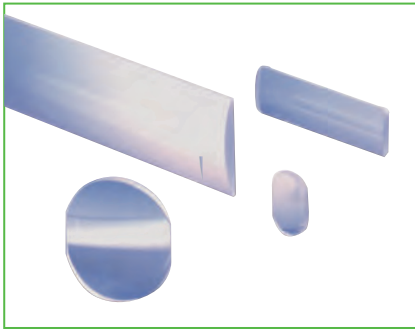Complex, optically active surfaces, which cannot be described by either spherical or aspheric equations, but are parameterized globally or locally by different mathematical models.

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
Cylindrical lenses

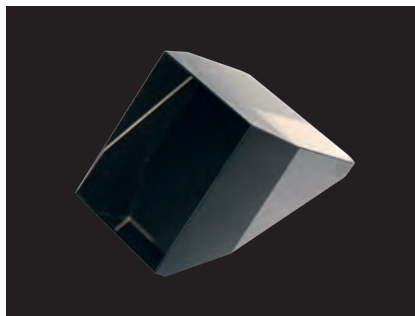- Rectangular, circular, and custom sizes
- Up to a maximum edge length of 0.4 m
- Custom coating design

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
Planar Optics (filters,prisms, tapered lenses)

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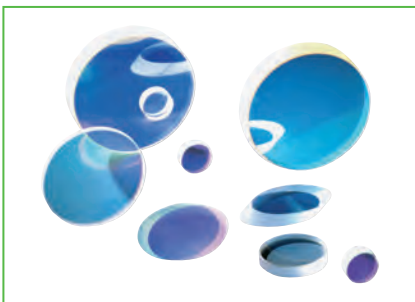
Coatings

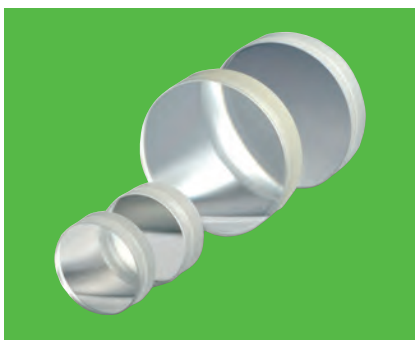- Antireflex coatings
 - From single coatings to
 - Ultra broadband coatings
- Band-pass filter
- Edge filter
- Mirror coatings fabricated from
 - Silver
 - Aluminum
 - Dielectric coatings

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Achromats

- All well-established cements
- Edge varnishing to reduce scattered light on request
- Diameter: Ø 3 mm up to Ø 80 mm

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CNC vertical integration:

- Small quantities (average lot size 20 pieces) in high quality
- Processing of all current glass materials incl. logging
- Glass technology laboratory for testing the optical glass values

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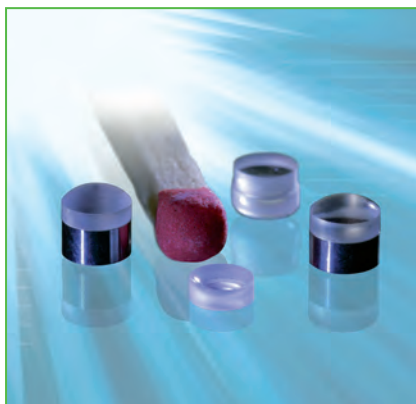
Micro Optics - Our Site in Asslar

The site in Asslar has been part of the Qioptiq Group since 2005, but its "optical competence" reaches much further back into the past. In 1952 it was founded as Neeb-Optik GmbH in Wetzlar, one of the German centers of optical industry. The company is specialized in the development and manufacture of micro optical systems and employs more than 80 skilled workers.

The production flow covers all work steps from grinding and polishing the lenses through coating up to micro assembly. The components produced are used not only in medical applications, in particular endoscopy, but also in many industrial fields, such as metrology and automation technology. The Asslar production site supplies both OEM (original equipment manufacturers) and business units within the Qioptiq Group, which further enhance these products by integrating them into complex systems.

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
Micro Optics, General

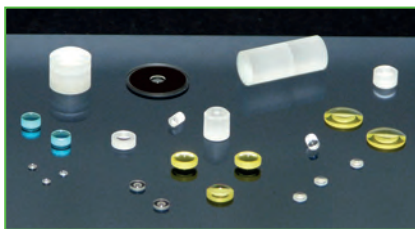- Grinding and polishing of spherical and plano optical components $\geq 0.3\text{mm}$ from all established optical lenses
- CNC and traditional production technologies
- Surface treatment suitable for laser quality
- Thin film coating of precision optics (AR, filters, mirrors and hydrophobic)
- Lithographic aperture stops and masks
- Optical fine cementing and alignment
- Assembling, testing and documentation
- Narrow tolerances for high end applications, e.g. 3D
- Wide range of Qioptiq test plates available in Zemax OpticStudio TPL
- Special test equipment

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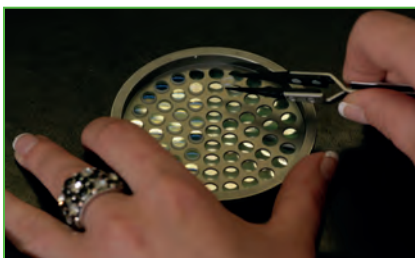
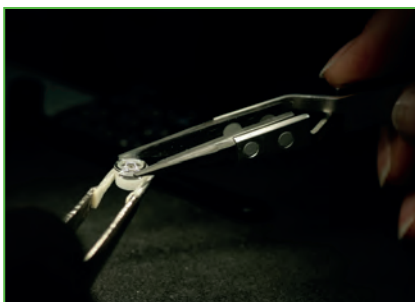
Customized Micro Optical Components

- Spherical lenses, doublets, triplets
- Rod lenses and compact objectives for rigid endoscopes
- Plano optics, e.g. cover glasses
- Optical glasses and fused silica
- Built-to-print according to customer drawing or specification
- From prototype to serial production

Spherical optics

- Lens diameter: 0.3 - 20 mm
- Possible surface quality: ≥ 0.5 fringes
- Possible irregularity: ≥ 0.2 fringes
- Centering error: $\geq 1'$
- Center thickness: up to ± 0.01 mm
- Diameter tolerance: ≥ 0.005 mm

Coating

- Single layer for all wavelengths
- Double layer (V-coating)
- Multi layer for UV to NIR
- Anti-reflective coating (e.g. MgF₂)
- Beamsplitter for specific wavelengths
- Back and front mirrors with Al and Ag
- Filters (e.g. IR-cut)

Cementing

- Diameter ≥ 0.5 mm (smaller diameters on request)
- Doublets – triplets – compact objectives – rod lens systems
- All typical UV and epoxy glues

Customized Opto-Mechanical Assemblies

- Development of optical systems according to customer-specific requirements
- Mechanical design
- Coating design
- Prototypes, pre-serial production and serial production
- Testing and documentation

Customized Micro Objectives

- Field of view: up to 170°
- Diameter: ≥ 0.8 mm
- Pinhole diameter: ≥ 0.1 mm ± 0.005 mm
- High precision field stop with chromium layer within the glued surface of the optical component, decentering ± 0.01 mm

Ingo Gretschel, Sr. Director Micro Optics and Factory Manager of our Asslar site

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Major tasks for small lenses

Originally, our expertise was mainly used in medical fields of application: endoscopy is one of the most advanced areas of optical miniaturization.

However, today our optical micro-lenses are needed in various industries and fields of application. Qioptiq is one of the leading providers of specialized micro-optics for industrial image processing and medical technology applications.

Our core business is the development, production and assembly of miniaturized optics. We produce spherical lenses, systems and optical assemblies to the highest level of precision and diameters from 0.3 mm to 15 mm. We process all optical glasses with classic fabrication methods.

Even if micro-optics continues to rely on conventional production methods, new technologies will expand applications of our products and improve their performance. Use of aspherical and GRIN lenses in our systems are examples of such opportunities.

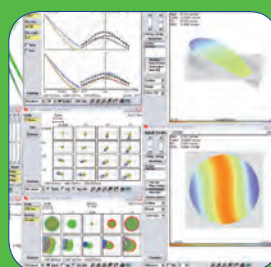
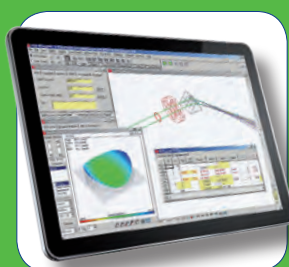
Optics Software

User-friendly Design Programs for Selection,
Analysis, and Optimization

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Our quality criteria:

- *Many years of experience in optical design and objective development*
- *Special design programs of WinLens solve your optical design challenges*
- *Easy-to-use selection software MachVis*
- *Continuous software development of both programs*

LINOS Optics Software

For nearly 20 years we have continuously developed, used and supplied WinLens - a user friendly and effective lens design program.

Our optical designers have decades of experience in modelling, analysis, optimisation and tolerancing - which feed into the ongoing development of this software suite. The simple interface provides an easy start for engineers - something which is just as important for training and teaching purposes.

Our LINOS MachVis software is designed to help you find the right lens for your machine vision application.

Starting with 4 simple parameters object size, working distance, sensor size, camera mount MachVis calculates the necessary optical parameters and then, from the extensive LINOS lens catalogue, selects all suitable lenses including matching accessories that meet the specifications.

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Special Features:

The WinLens Basic, PreDesigner, and MachVis programs are available for download free of charge. Databases of LINOS singlets, achromats, prisms and objectives are implemented.

Ideal Areas:

WinLens offers simulation, analysis and optimization of optical systems - both for training and teaching purposes as well as for professional use.

Inhalt

WinLens™

Softwarepaket WinLens™ 3D	488
Freeware	490
WinLens™ 3D	491
WinLens™ 3D Basic	491
Tolerancer	494
Glas Manager	496
Material Editor	498
PreDesigner	499
Lens Design Elements	502

MachVis

MachVis Software	500
MachVis Workflow	501

Content

WinLens™

<i>Software Package WinLens™ 3D</i>	<i>488</i>
<i>Freeware</i>	<i>490</i>
<i>WinLens™ 3D</i>	<i>491</i>
<i>WinLens™ 3D Basic</i>	<i>491</i>
<i>Tolerancer</i>	<i>494</i>
<i>Glass Manager</i>	<i>496</i>
<i>Material Editor</i>	<i>498</i>
<i>PreDesigner</i>	<i>499</i>
<i>Lens Design Elements</i>	<i>502</i>

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- WinLens™ 3D, Basic
- PreDesigner
- Lens Library

Software Paket WinLens 3D Software Paket WinLens 3D

Produktbezeichnung	Artikel-Nr.
Item Title	Part No.
Software package WinLens 3D Suite	G302020000

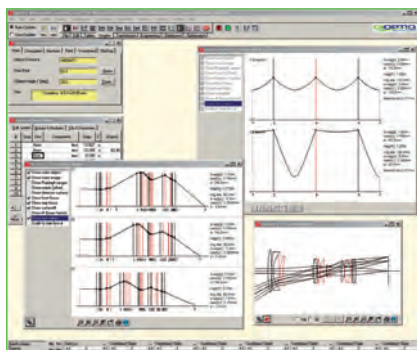
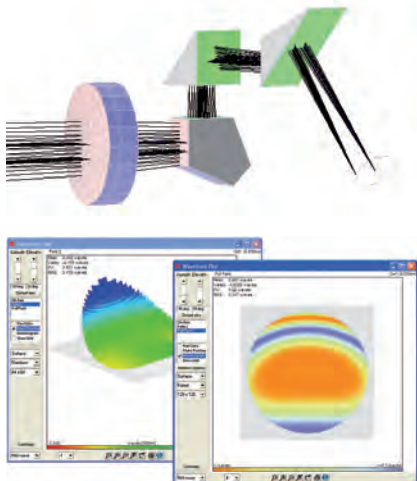
Software Package WinLens™ 3D

The software package can be used for PC's running Microsoft Windows™ and consists of four parts:

- WinLens™ 3D
- Tolerancer
- Glass Manager
- Material Editor


To use the software package WinLens™ 3D a license is required, which you receive by ordering this product. The following programs are available as freeware at www.winlens.de:

- WinLens™ 3D, Basic
- PreDesigner
- Lens Library

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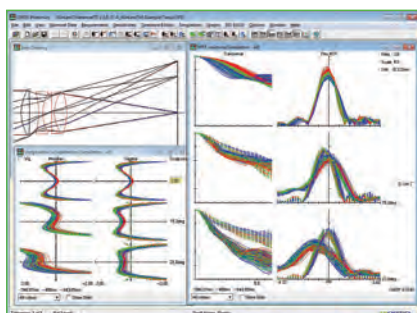
- Software for professional optical design
- Ideal tool for developing components and systems
- Meaningful charts and tables for analyzing optical systems
- Suitable both for our components (lenses, achromatic lenses, prisms, etc.) and for customer-specific components and systems
- Powerful optimization routine using relative defects
- Comprehensive engineering features
- Simulations of on-axis and folded systems

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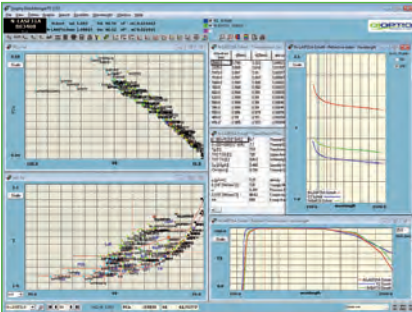
- For tolerancing the construction parameters of optical on-axis systems
- Statistical analysis of production tolerances
- Analyze the sensitivity of individual parameters
- Ideal tool for development and production of optical systems; can help reduce costs
- Create ISO 10110 drawings with highly flexible editor

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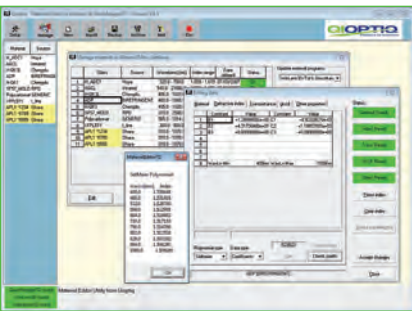- *Optical glass database of products from all leading suppliers*
- *Glass catalog data clearly laid out in tables and charts at standard or custom wavebands*
- *Sort and search for all types of glass*
- *Enter search criteria to select glass*
- *Vital tool for development and production of optical components and systems*

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- *Utility to create, edit & manage data for custom materials and melt derivatives*
- *Notes all key data in local files*
- *Simple to update the material databases in WinLens™ 3D, Tolerancer and Glass Manager*

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- *For more information, manuals, downloads and updates visit our website www.winlens.de*
- *For video tutorials visit www.opticalsoftware.net.*

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- Minimum PC configuration:*
- *PC running Windows 7, 8 or 10*
 - *Pentium Processor (300 MHz minimum), 64 MB RAM*
 - *Monitor resolution: ideally 1024x768 or better*
 - *Hard-drive installation space: 30 MB*

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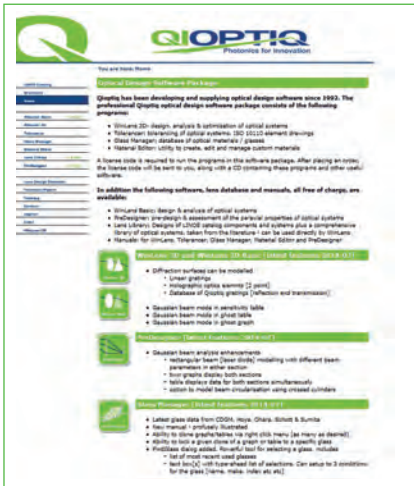
- *Fast: Run calculations right in lab*
- *Practical: Illustrated optical equations*
- *Mobile: For Android Mobiles and Tablets*

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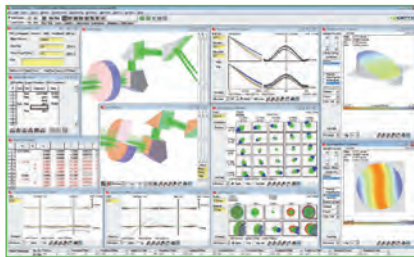
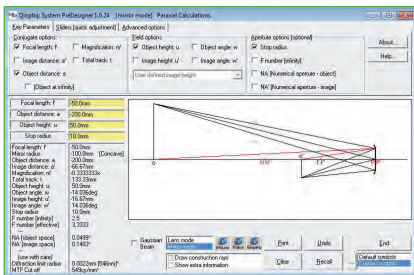
Freeware

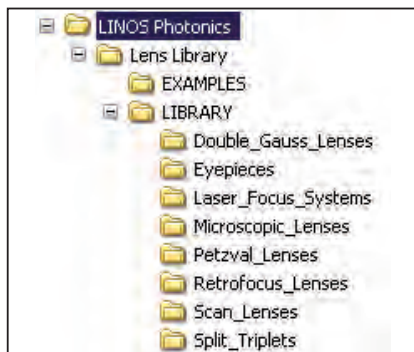
Minimum PC configuration:

- PC running Windows 7, 8 or 10
- Pentium Processor (300 MHz minimum), 64 MB RAM
- Monitor resolution: ideally 1024x768 or better
- Hard-drive installation space: 30 MB

In addition the following programs are available free of charge:

- Freeware version of the WinLens™ 3D optical design program
- For component and system development
- Informative charts and tables for analyzing optical systems
- Suitable both for our components (lenses, achromatic lenses, prisms, etc.) and for customer-specific components and systems
- Ideal educational tool for training in technical optics

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

- Free database of optical systems and special components (Qioptiq products) for use with WinLens™
- Additional comprehensive lens library with systems from specialized journals and patent literature
- Includes a selection of sample systems based on standard components

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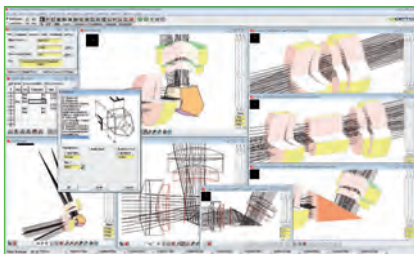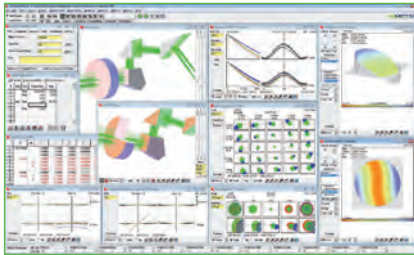* These functions are not available in the WinLens™ 3D Basic edition.

WinLens™ 3D WinLens™ 3D Basic

System

- Spherical and aspherical surfaces (rotationally symmetric, surface description also with standardized coefficients and Zernike coefficients)
- Cylindrical and toric surfaces
- Lenses, mirrors, blocks & prisms
- Surface coatings (Qioptiq coatings)
- Media between components (air, water, special media)
- Tilts and decenters for surfaces, components and groups
- Zoom systems (multiple configurations)
- Thin lens layout for pre-designs and starting systems (paraxial approach) *

Parameters

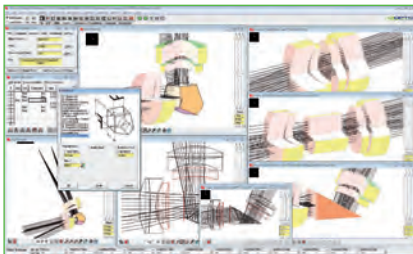
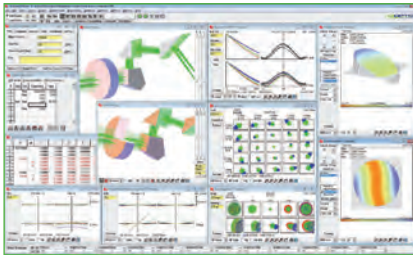
- Object and image positions
- Object and image sizes
- Magnification
- Aperture (numerical apertures, pupil radii, aperture radius, f-stop)
- Wavelengths
- Define media in object and image space (air, water, glass, special media)
- Radius of image curvature
- Defocussing

Analysis

- Analyses for any wavelength
- Lens drawing with beams (optional data output in DXF file)
- Basic optical data:
 - Paraxial data (focal point, cardinal points and pupil positions)
 - Paraxial ray tracing
- System data in table form
- Basic data on individual surfaces, elements and components
- Lens form analysis: ray heights, edge separations *
- Form analysis of aspherical surfaces *
- Lens weight
- Internal absorption loss
- Transmission
- Sensitivity analysis with variation tables for paraxial data and Seidel errors *
- Ray tracing
- Gaussian beam table for laser applications
- Image quality graphs:
 - Optical path differences (wave aberrations)
 - Transverse aberrations
 - Longitudinal aberrations
 - Chromatic aberration
 - Deviation from sine condition isoplanatism *
 - Field aberrations (curvature, astigmatism, distortion, color)
 - Spot diagrams
 - Gaussian graph
 - Geometric modulation transmission function (MTF):
 - MTF as a function of spatial frequency
 - MTF as a function of image size *
 - Through-focus MTF
 - Delano diagrams

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WinLens™ 3D WinLens™ 3D Basic

Analysis, continuation

- Relative illuminance *
- Footprint *
- Transmission functions (axial and for all chief rays)
- Seidel bar chart
- Ghost analysis *
- Graphical and tabular evaluation
- Interferogram and wavefront plot

Editing

- System data editor for components
- System data editing in surface tables
- Unlimited numbers of components and configurations (zooms)
- Linking of components (pick-ups)
- Tool fitting for existing Qioptiq tools
- Sliders and scaling of elements and groups
- System parameters (conjugates, aperture, field, wavelengths)
- Wavelength editor (for selecting standard and custom wavelengths)
- Selection of alternative glass types
- Data on real glass melts

Databases

- Qioptiq components (singlets, achromatic lenses, prisms, wedge plates and more)
- Choice of search criteria includes:
 - Lens type
 - Unmounted, mounted, Microbench mounted
 - Focal length range
 - Aperture range
- Material (optical lenses, special materials):
 - Databases of glasses from all leading suppliers
 - Interactive glass selection with customizable glass diagrams
 - User-defined material database
- Coating (Qioptiq coatings for optical surfaces):
 - Broadband anti-reflective coatings
 - Narrowband anti-reflective coatings
 - Adaptable center wavelength
- Optical systems (separate database/Lens Library):
 - Beam expanders
 - Condenser systems
 - Microprojection lenses
 - Large-aperture laser lenses
 - Microscope lenses, eyepieces
 - Comprehensive system library based on specialized and patent literature (separate download)

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Optimization *

- *Efficient optimization routines*
- *Choice of variable system parameters*
- *Choice of parameter limits*
- *Define merit function by relative defects*
- *Determine efficacy of variables*
- *Use special help routines to set up the merit function*
- *Replay individual optimization steps using the "Recorder" function*

Additional features

- *Sliders: Simple variation of system data to adjust individual components*
- *User-defined glass diagram*
- *Import/export data files to Code V, Oslo, Sigma & Zemax*
- *Export tables to Excel and HTML files*
- *Comprehensive representations, using:*
 - *2D or 3D lens drawing*
 - *Lens drawing/graphs can show data for selected zooms*
 - *Create multiple clones of drawings, graphs & tables*
 - *Individual scales for each clone*
 - *Freeze individual tables/graphs for system comparisons*
 - *Bookmark function for system comparisons*
- *Define modules and groups for subsystem analysis **

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* *These functions are not available in the WinLens™ 3D Basic edition.*

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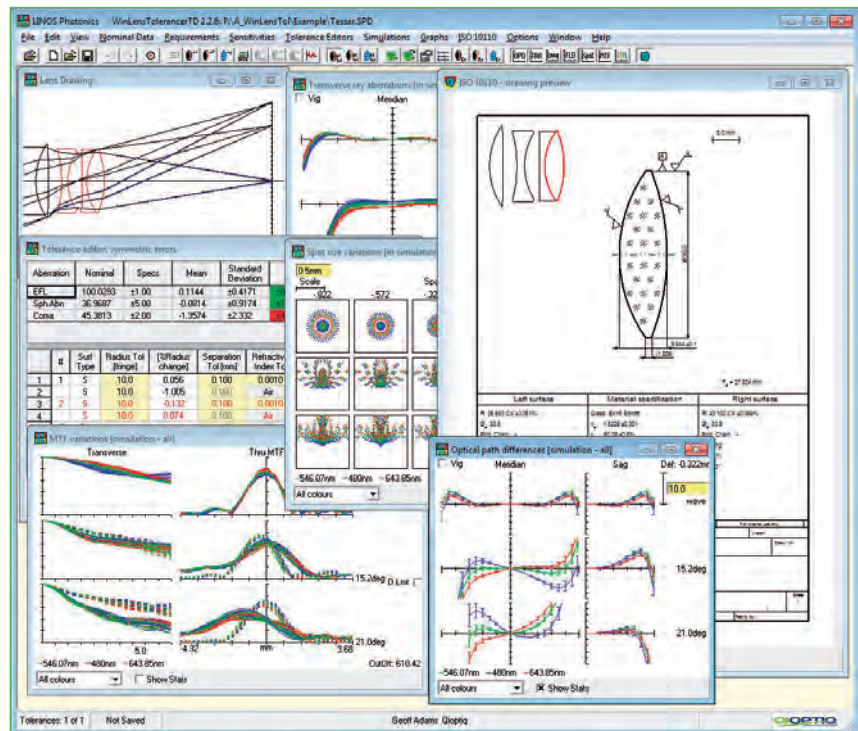
Our WinLens poster "Lens Design Elements" shows many helpful optical design illustrations, formulas, definitions and other information. The poster is shown on the last page of this chapter. Feel free to download it from www.winlens.de.

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Sensitivity analysis for symmetric, asymmetric and form errors

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Tolerance Editors

- *Simple definition of production tolerances in tables (centering errors, surface form deviations, etc.)*
- *Optional automatic tolerancing through a choice of default tolerances in several accuracy levels (normal, precise, extra precise)*
- *Simple definition of permissible performance variations in tables (paraxial data, Seidel aberrations, field aberrations, beam aberrations)*
- *Estimation of tolerance efficiency with respect to required maximum variations of paraxial data and aberrations*
- *Color-coded warning signals when permissible range of output data is exceeded*
- *Determination of tolerances using inverse sensitivity calculation*

Statistical tolerance analysis

- *Monte-Carlo analysis of optical systems with pre-defined production tolerances, for:*
 - *Determination of variation ranges*
 - *Compensation during simulation, for example through focus adjustment*
 - *Graphic display of aberrations and output charts during Monte-Carlo analysis*
 - *Graphic display of the standard deviation*
 - *Tabular comparison of nominal and Monte-Carlo simulated aberrations*
 - *Display performance of any system in the simulation*
 - *Display performance of a system with user defined errors*
- *ISO 10110 drawings*
 - *choice of english or german marking*
 - *customisable symbols*
 - *Output via dxf or bmp*

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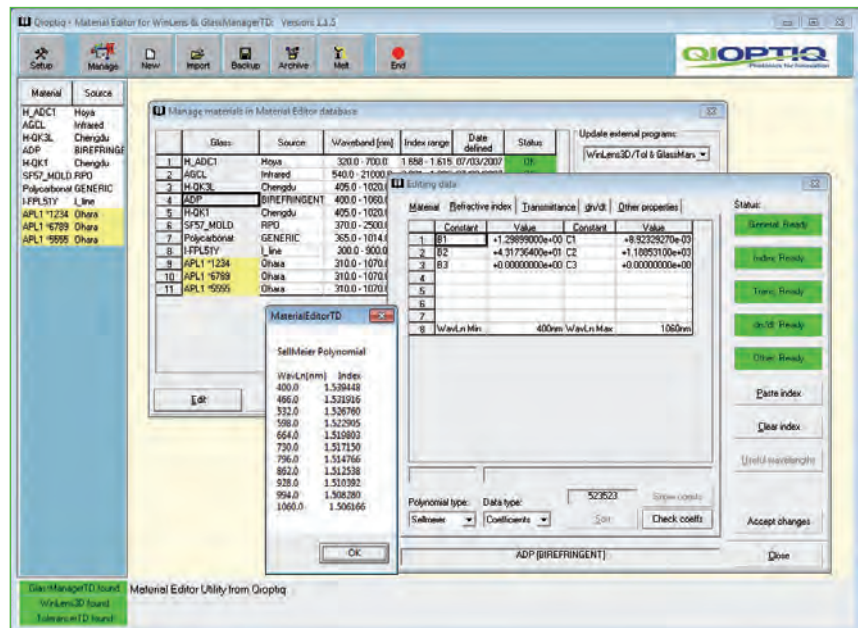
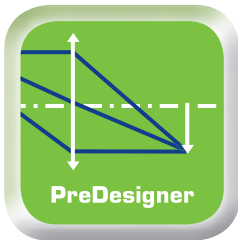
The mobile app for optical designers on the go! Convenient: Paraxial calculations via touchscreen

- Fast: Run calculations right in lab
- Practical: Illustrated optical equations
- Mobile: For Android Mobiles and Tablets

Download now for free in the Google Play Store !

MachVis Software

MachVis Software

- Video Tutorials finden Sie unter www.opticalsoftware.net.
- For video tutorials visit www.opticalsoftware.net.

Qioptiq Lens Name	FOCUS Device	Exit	Lens EFL	Total W.D.	Free W.D.	Lens to Camera	Flange F.D.	Flange F.D. Range
Theory			96.77mm	620.00mm	500.00mm		120.00mm	
1 Optem 0 24x 0 4xLL 0 6xCT	Absocal	0	Absocal	490.00mm			105.00mm	
2 Rogonar-S 90	Modular Fo.	1	90.10mm	578.44mm	446.12mm	114.82mm	101.52mm	97.7-122.7mm
3 Apo-Rodagon N 90	Modular Fo.	1	90.10mm	574.04mm	436.32mm	120.22mm	106.52mm	97.7-122.7mm
4 Rogonar-S 90	Smart Focus	2	90.10mm	578.44mm	446.12mm	114.82mm	101.52mm	93.2-105.6mm
5 Apo-Rodagon N 90	Smart Focus	2	90.10mm	574.04mm	436.32mm	120.22mm	106.52mm	105.2-117.6mm
6 Apo-Rodagon N 90	Focus Adap.	7	90.10mm	574.04mm	436.32mm	120.22mm	106.52mm	104.5-109.5mm
7 Apo-Rodagon N 105	Modular Fo.	2	105.20mm	670.68mm	510.23mm	142.95mm	124.45mm	122.2-147.2mm
8 Apo-Rodagon N 105 full metal	Modular Fo.	2	105.20mm	670.68mm	510.23mm	142.95mm	124.45mm	122.2-147.2mm
9 Apo-Rodagon N 105	Smart Focus	3	105.20mm	670.68mm	510.23mm	142.95mm	124.45mm	117.2-129.6mm
10 Apo-Rodagon N 105 full metal	Smart Focus	3	105.20mm	670.68mm	510.23mm	142.95mm	124.45mm	117.2-129.6mm
11 Rodagon 105	Modular Fo.	1	106.40mm	660.77mm	523.03mm	140.24mm	127.04mm	125.2-150.2mm
12 Rodagon 105 azimuth	Modular Fo.	1	106.40mm	660.77mm	523.03mm	140.24mm	127.04mm	125.2-150.2mm
13 Rodagon 105 full metal	Modular Fo.	1	106.40mm	660.77mm	523.03mm	140.24mm	127.04mm	125.2-150.2mm
14 Rodagon 105	Smart Focus	3	106.40mm	660.77mm	523.03mm	140.24mm	127.04mm	117.2-129.6mm
15 Rodagon 105 azimuth	Smart Focus	3	106.40mm	660.77mm	523.03mm	140.24mm	127.04mm	117.2-129.6mm
16 Rodagon 105 full metal	Smart Focus	3	106.40mm	660.77mm	523.03mm	140.24mm	127.04mm	117.2-129.6mm
17 Rodagon 105	Focus Adap.	0	106.40mm	660.77mm	523.03mm	140.24mm	127.04mm	126.5-131.5mm
18 Rodagon 105 azimuth	Focus Adap.	0	106.40mm	660.77mm	523.03mm	140.24mm	127.04mm	126.5-131.5mm
19 Rodagon 105 full metal	Focus Adap.	0	106.40mm	660.77mm	523.03mm	140.24mm	127.04mm	126.5-131.5mm

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Starting with 4 simple parameters:

- Object size
- Working distance
- Sensor size
- Camera mount

MachVis calculates the necessary optical parameters and selects all suitable lenses from the extensive LINOS machine vision portfolio that meet the specifications.

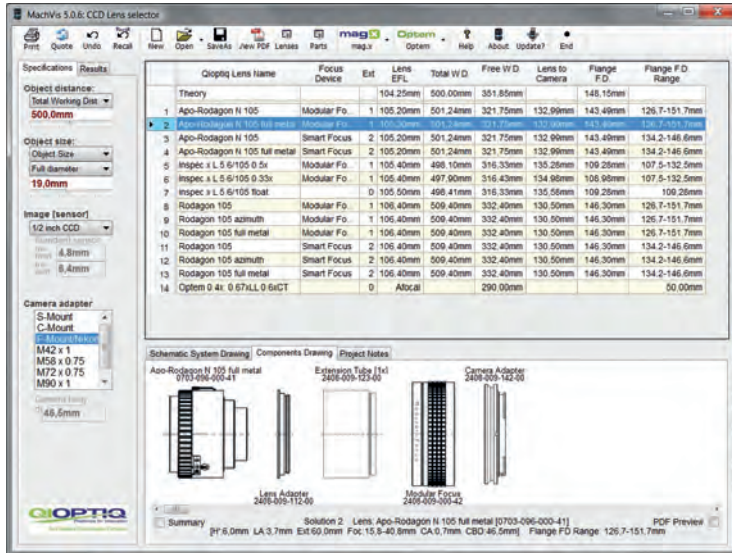
If mechanical accessories (e.g. focussing unit, extension tubes etc.) are needed, MachVis creates a printable, itemised list of required parts. MachVis also generates a schematic drawing of the lens and accessories in the order of assembly.

Specifications, along with your own project notes, may be saved to disk for later recall. Datasheets (PDF) for each lens may be viewed for a more detailed analysis.

- Paraxial geometric optic
- Automatic selection of suitable systems
- Lens database
- Mechanics database
- Data sheets and 3-D models
- Configurator for mag.x (high-resolution wide-field microscope)
- Configurator for Optem Fusion

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
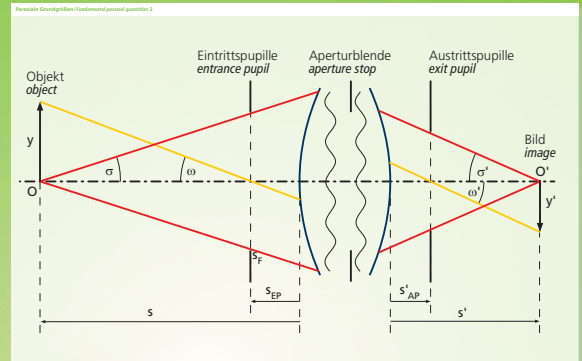
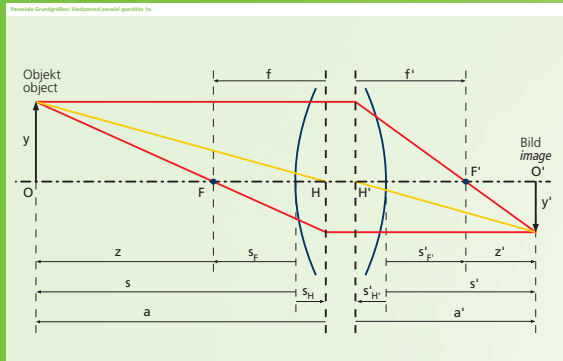
For download and further information, please go to www.qioptiq.com/request-machvis-software.html.

Download now for free!

LENS DESIGN ELEMENTS

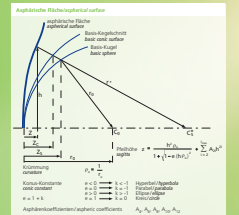Grid of ray diagrams for various lens types: plano-convex, biconvex, meniscus, aspherical, cylindrical, etc.

Grid of ray diagrams for more complex lens systems like microscope, telescope, and camera lenses.

Table with optical formulas: Snell's law, lens maker's equation, magnification, etc.

Table with paraxial optics formulas: Gaussian beam parameters, ABCD matrices, etc.

Table with aperture formulas: NA, f-number, etc.

Table with dispersion formulas: Sellmeier equation, Abbe number, etc.

Table with field geometry formulas: field of view, etc.

Table with diffraction limit formulas: Rayleigh, Sparrow, etc.

Table with Gaussian beam formulas: waist, divergence, etc.

Table with aberration formulas: spherical, coma, astigmatism, etc.

Table with wavelength range data: IR, VIS, UV, EUV ranges.

Table with symbol key: a, a', n, f, etc. symbols used in diagrams.

Table with image format data: Type, Format, Diagonal, etc. for various camera formats.

Table with wellenlänge / wavelength data: Laser types and their wavelengths.

Table with wellenlänge / wavelength data: Laser types and their wavelengths.

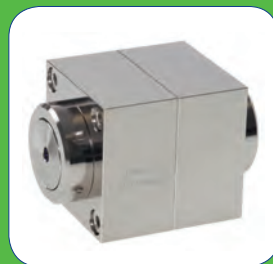
Electro Optics

High-performance Faraday isolators,
Pockels cells and modulators

Faraday Isolators

Laser Modulators and Pockels Cells

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Our quality criteria:

- Isolation ≥ 30 dB (one-stage) or ≥ 60 dB (two-stage)
- Transmission ≥ 90 % (one-stage) or ≥ 80 % (two-stage)
- All models can be used over wide wavelength ranges

LINOS Faraday Isolators

We have Faraday isolators for all wavelengths in the range from 390 to 1310 nm, as well as 1550 nm. Isolators for other wavelengths can be offered upon request. Many isolators can be adjusted over a wide spectral range; tunable frequency models can even be set for an interval of several hundred nanometers. At the same time, LINOS Faraday isolators are distinguished by high performance combined with the greatest possible transmission. With more than 60 dB, our two-stage isolators offer the best isolation available on the market.

The consistently-high Qioptiq quality is assured by a combination of our many years of experience, an sophisticated design, modern engineering with computer simulations and sophisticated processing. The result is the outstanding value that distinguishes all our products – value you can count on!

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Special Features:

Our isolators are available in many different versions for a large range of wavelengths.

Ideal Areas:

Protecting lasers from damage or instability; Decoupling oscillators and amplification systems; Injection locking panels, and more.

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Content

<i>Characteristics and Technical Features</i>	510
<i>Overview Faraday Isolators</i>	511
<i>Single Stage Faraday Isolators, Technical Overview</i>	512

Faraday Isolators

<i>Isolators with 2 mm Aperture, SV Series</i>	513
<i>Isolators with 3.5 and 5 mm Aperture, SC Series</i>	514
<i>Low outgassing Isolators with 3.5 mm aperture, LO Series</i>	516
<i>Isolators with 5 mm Aperture, SV Series</i>	517
<i>Isolator with 8 mm Aperture, SI Series</i>	518
<i>Special Isolators with 5 mm Aperture, SV Series</i>	519
<i>Isolators with Magneto-Optical Crystal Film</i>	520
<i>Isolators with a Broad Tuning Range</i>	521
<i>Tunable Isolators with 5 mm Aperture, SV/SI Series</i>	522
<i>Tunable Isolator with 8 mm Aperture, SI Series</i>	523
<i>Two Stage Faraday Isolators, Technical Overview</i>	525
<i>Two-Stage Faraday Isolators, tunable, TVC/TIC Series</i>	526
<i>Two-Stage Faraday Isolators, tunable, DLI Series</i>	527

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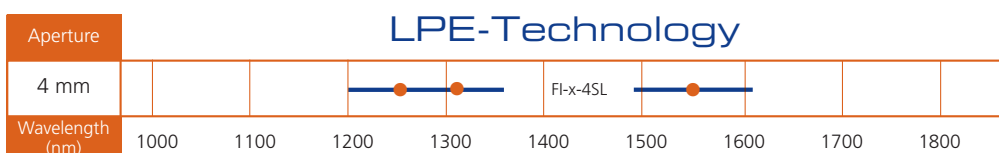
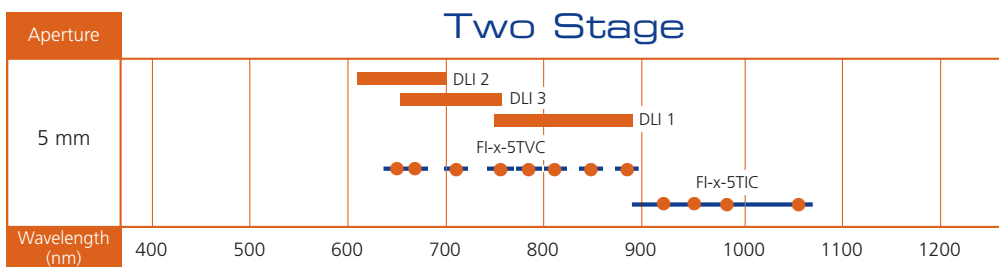
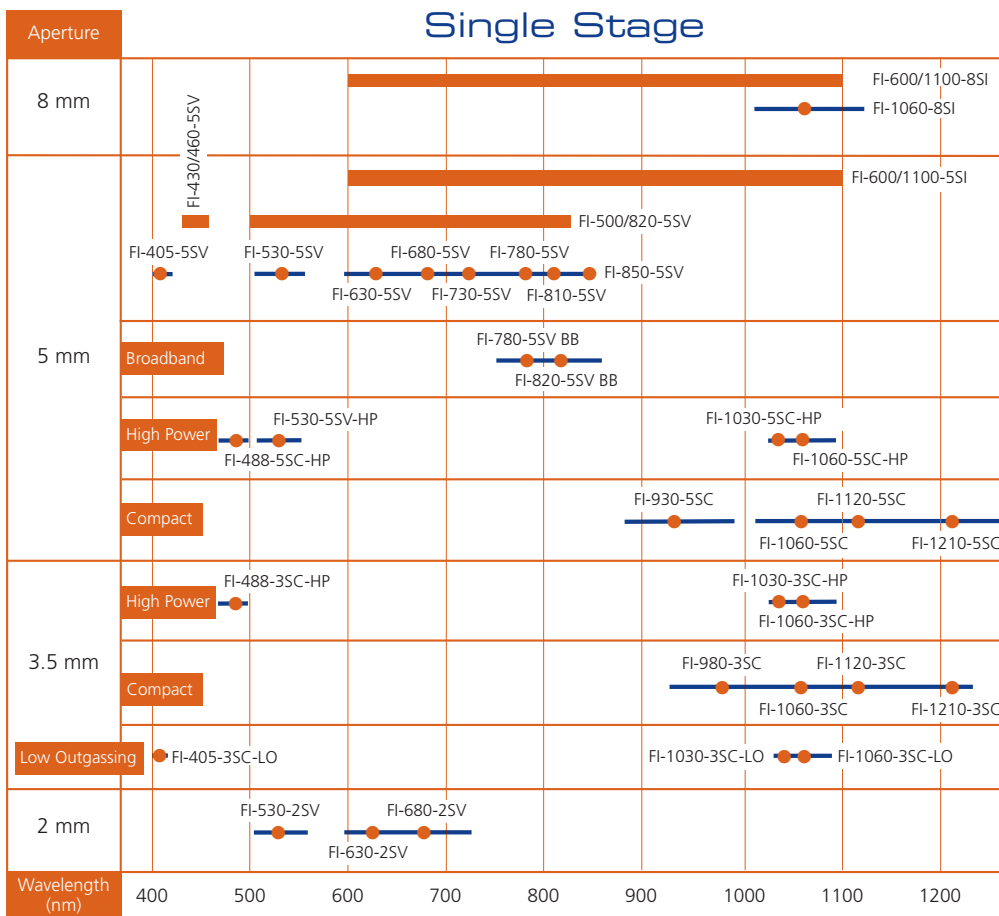
Characteristics and Technical Features

Principle of Operation

Faraday isolators are optical components which allow light to travel in only one direction. Their mode of operation is based on the non-linear Faraday effect (magneto rotation). In principle, the function of an optical isolator is analogue to that of an electrical diode.

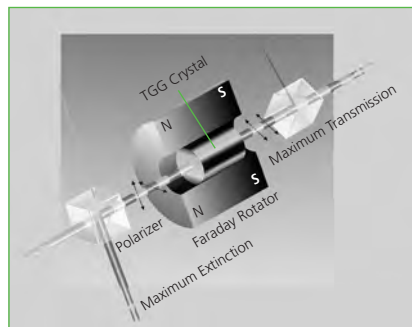
Faraday isolators are composed of three elements:

- Entrance Polarizer
- Faraday Rotator
- Exit Polarizer

Thin film polarizers are commonly used as entrance and exit polarizers, typically in form of a special polarizing beam splitter cube. These polarizers have an extremely high extinction ratio and are designed for use with high power lasers. The polarizer entrance and exit surfaces are coated with an antireflective coating for the specified wavelength range. The key element of the Faraday isolator is the Faraday rotator. The rotator consists of a strong permanent magnet containing a crystal with a high Verdet constant.

Light of any polarization entering the entrance polarizer exits as horizontally or vertically linearly polarized light. Since laser light is usually linearly polarized, one can match the orientation of the entrance polarizer and the laser by simply rotating the isolator. Light then passes through the Faraday rotator. For most wavelengths the crystal is a Terbium Gallium Garnet (TGG) crystal which is placed in a strong homogeneous magnetic field. Crystal length and magnetic field strength are adjusted so that the light polarization is rotated by 45° on exiting the crystal. In the figure above the light is rotated counter clockwise when viewed in the north/south direction of the magnetic field ($\pm 45^\circ$) and the exit polarizer is also oriented at $\pm 45^\circ$, so that the maximum beam intensity is transmitted.

If light of any polarization, but with a reversed direction of propagation, meets the exit polarizer, it leaves at $\pm 45^\circ$, passes through the Faraday rotator and is again rotated by $\pm 45^\circ$. The non-reciprocal nature of the Faraday effect results in the direction of rotation once again being counter clockwise as viewed in the north/south direction of the magnetic field. Upon leaving the Faraday rotator, the polarization has gone through two $\pm 45^\circ$ rotations resulting in a total rotation of $\pm 90^\circ$. In this polarization direction the light is deflected laterally by the entrance polarizer.

Increased Isolation

The maximum isolation of the Faraday isolator is limited by inhomogeneities of the TGG crystal and the magnetic field. However, it is possible to square the extinction ratio by placing two isolators in series and by arranging the polarity of the two magnets to be opposite to each other. This way the polarization direction of the transmitted light remains unchanged in the transmission direction and the effect of both magnetic fields is enhanced. This arrangement also leads to a more compact isolator.

Literature:

1. Bergmann, Schäfer
Lehrbuch der Experimentalphysik,
Band 3, Optik de Gruyter, Berlin 1993
2. R. Wynands et. al.
A compact tunable 60-dB Faraday
optical isolator for the near infrared
Rev. Sci. Instrum. 63, 5586-5590 (1992)
3. M. Lenzner et. al.
Sub-20-fs, kilohertz-repetition-rate
Ti:sapphire amplifier, Optics Letters,
Vol. 20, No. 12, 1397-1399

High Isolation

The properties of the Faraday isolator are determined by the quality of the optical elements and the uniformity of the magnetic field. The entrance and exit polarizers exhibit a very high extinction ratio, so that the isolation is mainly limited by inhomogeneities in the crystal material. Specially selected crystal materials with a high Verdet constant combined with permanent magnets with a high remanence enable us to use shorter crystals and obtain an isolation > 30 dB.

The radiation blocked by the entrance and exit polarizers is not absorbed internally, but is deflected by 90° with respect to the beam direction. This ensures a stable thermal operation even at higher laser power levels. The blocked radiation can be used for other applications. All optical surfaces are slightly tilted relative to the beam axis.

Low Insertion Loss

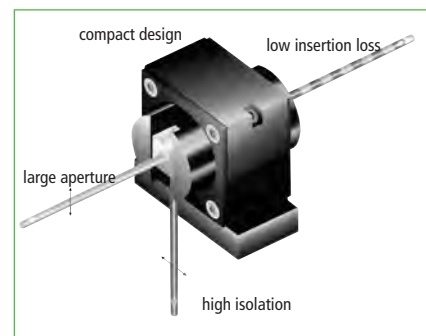
The high transmission, typically $> 90\%$, is achieved by using low absorbing materials and antireflective coatings with low residual reflectivity on all entrance and exit surfaces.

Large Aperture, Compact Design

All optical elements have been aligned to eliminate beam shading and allow for easy adjustment. Focusing is not necessary.

The compact design is achieved by using rare earth magnets with the highest remanent magnetism and TGG crystal material with a high Verdet constant. The isolator is suitable for divergent beams or in setups with limited space. A minimal optical path length in the isolator results in the lowest possible influence on the image.

Three sides of the entrance and exit polarizers are usable and readily accessible for easy cleaning. The degree of isolation can be adjusted in a wide range.

Mounting Flexibility

The isolators can be mounted directly via threaded holes in the housing or via additional base plates or angle brackets.

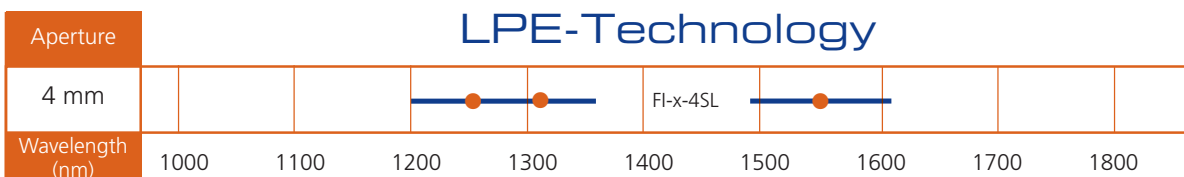
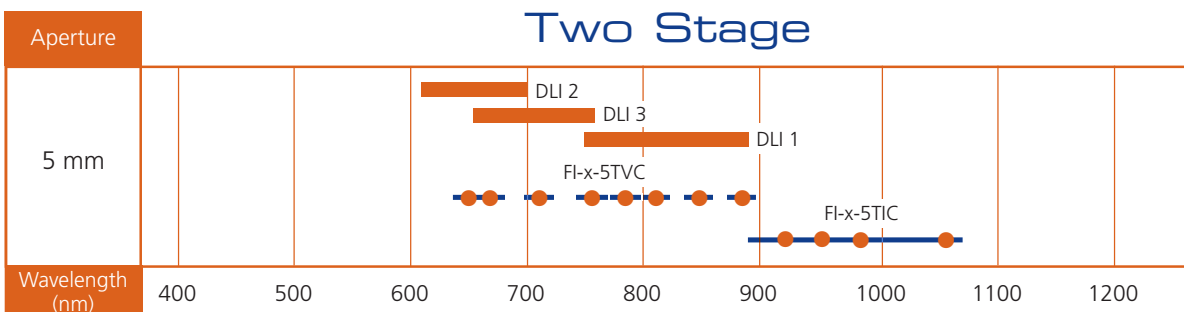
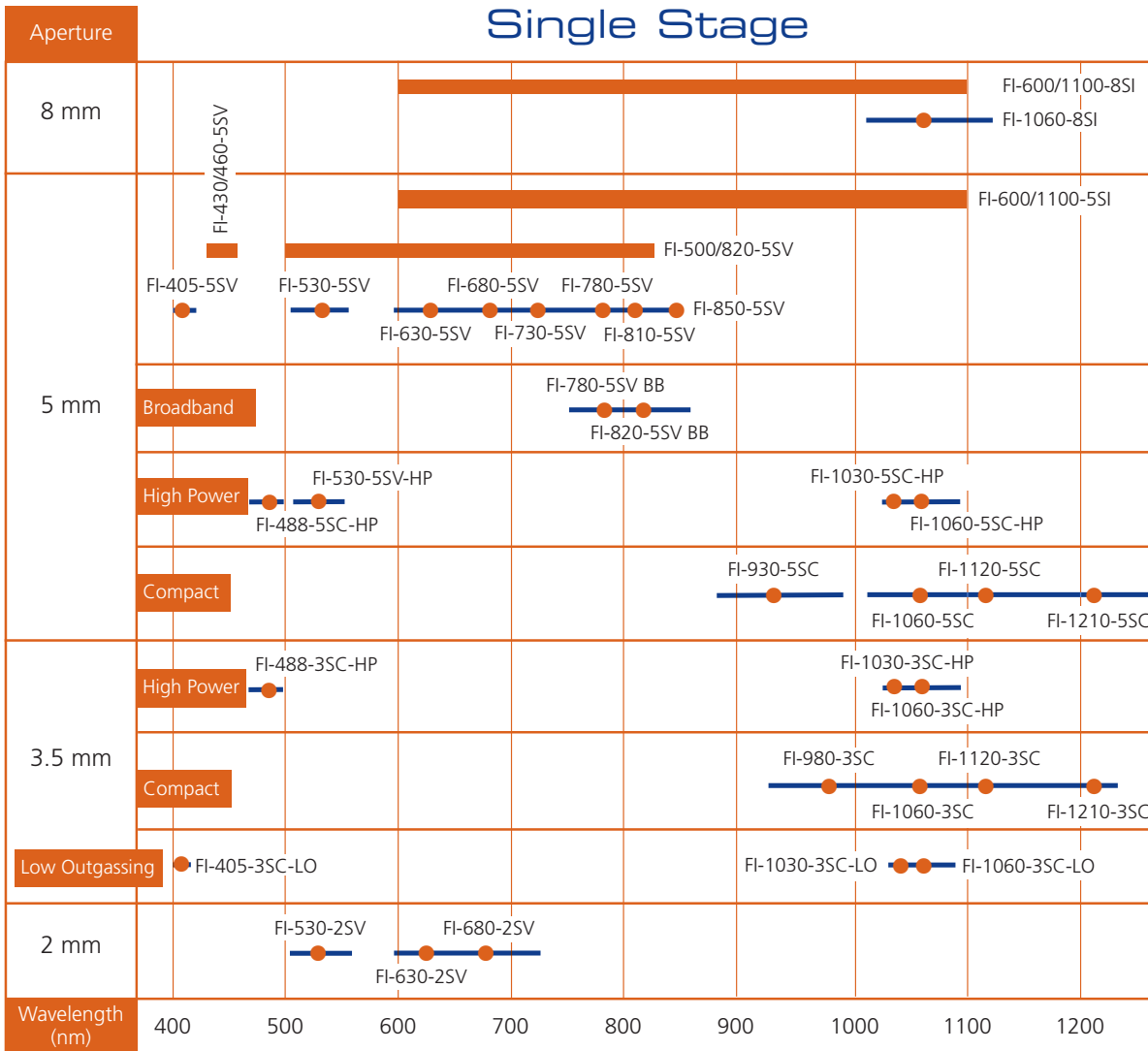
Applications

The ongoing development and refinement of laser technology have created a need for optical components that shield the laser resonator from back reflections. Faraday isolators provide an efficient method of suppressing instabilities and intensity fluctuations in laser devices.

Typical applications are:

- Protection of the resonator in solid state and gas lasers from back reflections
- Prevention of parasitic oscillation in multistage solid state amplifiers
- Protection of diode lasers against back scatter and extraneous light

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Single Stage Faraday Isolators, Technical Overview

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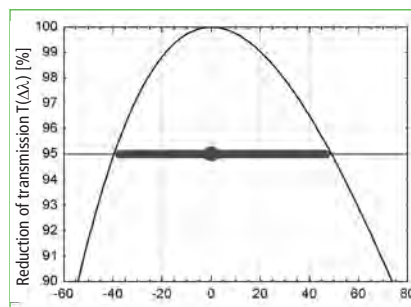
Tuning of design wavelength $\Delta\lambda = \lambda - \lambda_0$

Introduction

The compact Faraday isolators in this chapter use a single stage rotator. The length is kept to a minimum with the use of powerful permanent magnets in an optimized geometry. A 360° rotation of the exit polarizer provides a maximum extinction over a certain range around the central wavelength. The entrance and exit polarizers are polarizing beam splitter cubes. The blocked radiation is deflected by 90° and is readily available for other applications. At 30 dB, the specification of the isolator is sufficient for most standard applications. For specialized applications, selected isolators with an extinction up to 45 dB are available.

An even higher extinction is provided by the two stage isolator series.

Wavelength tuning

The Verdet constant of the TGG crystal is dependent upon wavelength and temperature. In order to compensate for different temperatures or different wavelengths, it is possible to tune the isolator in order to achieve maximum extinction. Tuning the isolator is accomplished by rotating the holder of the exit polarizer with respect to an engraved angular scale.

The graph shows the typical reduction factor of the transmission $T(\Delta\lambda)$ that is due to the tuning of the isolator to a wavelength λ that is different from the design wavelength λ_0 . The bar has a length that covers the wavelength range for which $0.95 < T(\Delta\lambda) < 1$. The bullet indicates the design wavelength λ_0 . The overall transmission of a Faraday isolator is equal to $T_t = T_0 \times T(\Delta\lambda)$, where T_0 is a factor that represents the intrinsic transmittance of the isolator. At the design wavelength the overall transmission of the Faraday isolator is $T_0 > 90\%$.

Broadband option (Special isolators)

On Broadband (BB) models the isolation is improved over a broadband spectrum by compensating the rotational dispersion of the TGG. This renders the device usable over a wavelength range of ± 50 nm without additional adjustment. The isolators can be mounted on rods, cylindrical mounts or by using the assembly surfaces so that the laser polarization can be oriented horizontally or vertically.

Applications

The following single stage Faraday isolators are suitable for all kind of lasers, especially:

- Ar+ and Kr+ lasers
- other Ion lasers
- HeNe lasers
- other Gas lasers
- Dye lasers
- Diode lasers
- Ti: Sapphire lasers
- Cr:LiCAF lasers
- Short Pulse lasers
- Mode-synchronized lasers
- Alexandrite lasers

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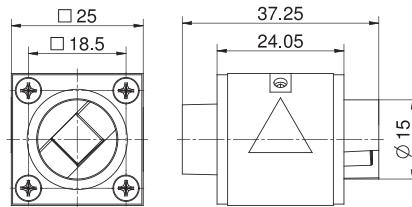
Isolators with 2 mm Aperture, SV Series

FI-x-2SV

- Isolation better than 30 dB / typically 38-42 dB over the entire wavelength range
- Custom isolation values on request
- TGG crystal
- Output polarizer, 360° rotation, engraved tuning scale
- Access to blocked beam
- Aperture Ø 2 mm
- Dimensions 25x25x37 mm³
- Mounting 2SV-version: via two M3 threaded holes at the bottom side, 20 mm separation

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Isolators with 3.5 and 5 mm Aperture, SC Series

FI-x-3SC / FI-x-5SC

- Extremely compact design
- Isolation better than 30 dB, typically 38-42 dB over the entire wavelength range
- TGG crystal
- Access to blocked beam
- Brewster polarizers available (BP)
- HP versions include optically contacted polarizers
- Integrated waveplate optional
- Mounting: via four M2 threaded holes on the bottom side and backside

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

Isolators series SC feature small and compact design with IR wavelength range.

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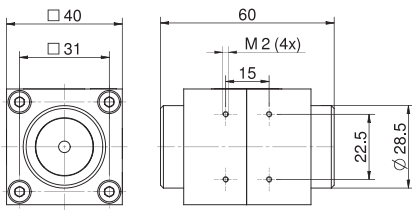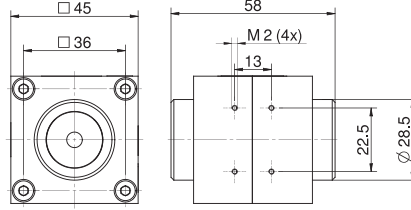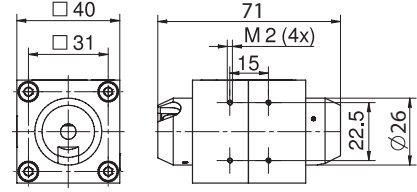
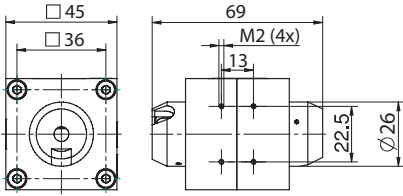
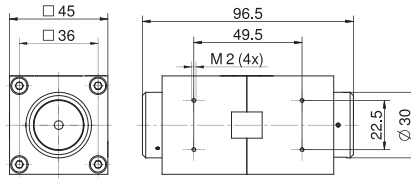
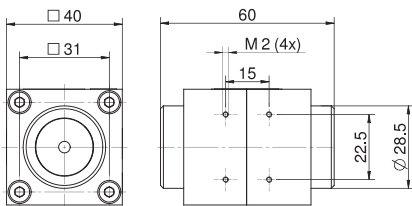
Low outgassing Isolators with 3.5 mm aperture, LO Series

FI-x-3SC LO

Our extensive experience and innovations in optical solutions for semiconductor manufacturing have led to the first ever Low Outgassing Faraday Isolators. These Isolators are ideally suited for integration in encapsulated and high-power laser systems in the wavelength range from the violet to the NIR. The unique design emits 25-times less volatile organic compounds (VOCs) than conventionally produced isolators.

Optically contacted, fused-silica input and output polarizers ensure high damage threshold while delivering highest isolation of 33 dB and outstanding transmittance of 95% in the NIR and 86% in the violet.

They feature a 3.5 mm clear aperture and a patented magnet system, which enables an exceptionally compact design footprint.

- Emitting 25-times less volatile organic compounds (VOC) than conventionally produced isolators
- Sophisticated, low outgassing design and assembly
- Isolation ≥ 33 dB, typically 38-42 dB over the entire wavelength range
- Extremely compact package
- Optically contacted polarizers for high damage threshold and long lasting quality
- Access to blocked beam
- Stable and flexible mounting via four M2 threaded holes on the bottom side and backside
- Custom wavelengths and specifications upon request
- Integrated waveplate optional

Ausgasarme Isolatoren mit 3.5 mm Apertur, LO-Serie Low outgassing Isolators with 3.5 mm aperture, LO Series

Produktbezeichnung	Isolation garantiert/typisch (dB)	Transmission bei den Designwellenlängen (%)	Transmission an den Eckwellenlängen (%)	Einstellbereich (typisch) (nm)	Apertur \emptyset (mm)	Abmessungen (mm)	Zerstörschwelle @ 1.2 ps Pulse (J/cm ²)	Zerstörschwelle @ 8 ns Pulse (J/cm ²)	Part No.
Item Title	Isolation, guaranteed/typical (dB)	Transmission at design wavelength (%)	Transmission at boundary wavelength (%)	Tuning range typical (nm)	Aperture \emptyset (mm)	Dimensions (mm)	Damage threshold @ 1.2 ps pulses (J/cm ²)	Damage threshold @ 8 ns pulses (J/cm ²)	Part No.
FI-405-3SC LO	> 33/38-42	≥ 86	≥ 86	400 - 410	3.5	40x40x60	0.2 @ 405nm	2.6 @ 488 nm	845110900026
FI-1030-3SC LO	> 33/38-42	≥ 95	≥ 90	1025 - 1064	3.5	40x40x60	0.6 @ 1030 nm	8.2 @ 1030 nm	845110100230
FI-1060-3SC LO	> 33/38-42	≥ 95	≥ 90	1030 - 1090	3.5	40x40x60	0.6 @ 1060 nm	8.5 @ 1060 nm	845110100231

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

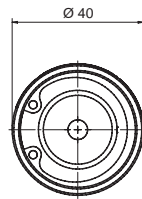
Isolators with 5 mm Aperture, SV Series

FI-x-5SV

- Isolation better than 30 dB, typically 38-42 dB over the entire wavelength range
- Custom isolation values on request
- TGG crystal
- Output polarizer, 360° rotation, engraved tuning scale
- Access to blocked beam
- Aperture Ø 5 mm
- Optional version with optically contacted polarizers on request, isolation better than 30 dB, typically better than 38-42 dB
- FI-405-5SV optically contacted polarizers as standard
- For upgrading to broadband-version refer to chapter Special Isolators
- Mounting: via two M4 threaded holes on the bottom side and the back side, 30 mm separation (5SV-version), or via base plate
- Base plate (50x30x9.5 mm³) included

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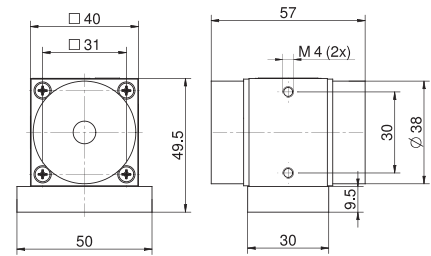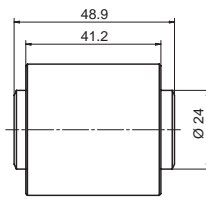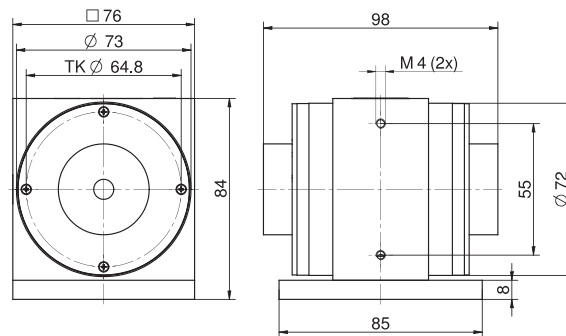
Isolator with 8 mm Aperture, SI Series

FI-1060-8SI


- Isolation better than 30 dB, typically 38-42 dB over the entire wavelength range
- Custom isolation values on request
- TGG crystal
- Output polarizer, 360° rotation, engraved tuning scale
- Access to blocked beam
- Optional version with Brewster plate polarizers (BP) on request, isolation better than 30 dB
- Optional version with optically contacted polarizer on request, isolation better than 30 dB, typically better than 38-42 dB
- Mounting: via two M4 threaded holes on the bottom side and the back side, 55 mm separation, or via base plate
- Dimensions 76x76x98 mm³
- Base plate (85x76x8 mm³) included

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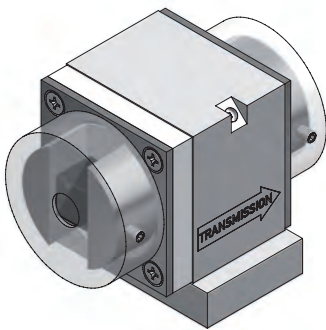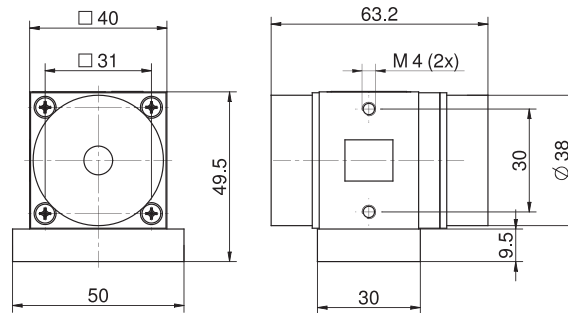
Special Isolators with 5 mm Aperture, SV Series

FI-x-5SV-MB / FI-x-5SV-BB

- Stable isolation better than 30 dB over the entire wavelength range without further adjustment
- Custom isolation values on request
- TGG crystal
- Output polarizer, 360° rotation, engraved tuning scale
- Access to blocked beam
- For multiline lasers or other broadband lasers such as fs-laser systems
- Mounting: via two M4 threaded holes at the bottom side and the back side; 30 mm separation, or via base plate
- Dimensions 40x40x63.2 mm³
- Base plate (50x30x9.5 mm³) included

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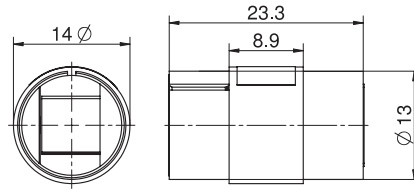
Isolators with Magneto-Optical Crystal Film

FI-x-4SL

- Extremely small size
 - Isolation better than 35 dB
 - Faraday Material: magneto-optical crystal film in saturation
 - Output polarizer, 360° rotation
 - Access to blocked beam
- Dimensions Ø 14x23.3 mm
 - Max. cw power: 8 W

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Isolators with a Broad Tuning Range

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Introduction

The function of the tunable isolators in the following chapter is based on a single stage Faraday isolator. Precision mechanics allow a continuous adjustment of the interaction between the magnetic field and the TGG crystal without moving any optical components.

It is possible to set the rotation angle to any value between 0° to 45° within the wavelength range in order to study the effects of varying degrees of feedback. Easy access to the blocked beam is provided by polarizing beam splitter cubes, which deflect the blocked beam by 90° .

Precision mechanics allow the exact reproduction of adjustments previously established. And with the addition of an optional micrometer display, an angular resolution in the arc minute range is achievable. The incorporation of very powerful magnets ensures a compact and efficient design.

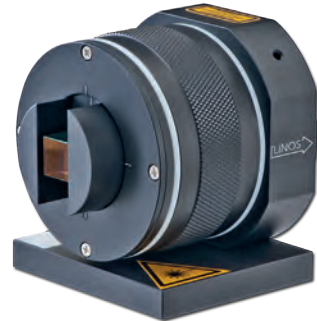
Operation

The isolator can be mounted on rods, cylindrical mounts or by using the assembly surfaces so that the laser polarization can be oriented horizontally or vertically. The entry and exit polarizers can be easily cleaned by removing the safety caps.

Applications

These isolators are suitable for all lasers operating in the 430-460 nm, 500-820 nm, respectively in the 600-1100 nm wavelength range especially for:

- Ar+ and Kr+ lasers
- other Ion lasers
- HeNe lasers
- Other gas lasers
- Diode lasers
- Nd:YAG lasers
- Ti: Sapphire lasers
- Cr:LiCAF lasers
- Dye lasers
- Alexandrite lasers
- Mode-locked lasers
- Short Pulse lasers

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
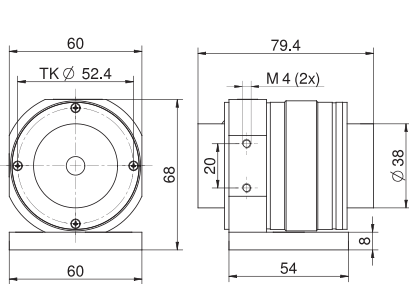
Tunable Isolators with 5 mm Aperture, SV/SI Series

FI-x/y-5SV / FI-x/y-5SI

- Continuous adjustment for wavelength without movement of optical components
- Tunable with maximum transmission and isolation over the complete wavelength range
- Isolation better than 30 dB, typically 38-42 dB over the entire wavelength range
- Custom isolation values on request
- TGG crystal
- Access to blocked beam
- Aperture Ø 5 mm
- Mounting: via two M4 threaded holes on the bottom side and the back side, 20 mm separation (5SV-version), 55 mm separation (5SI-version), or via angle bracket (5SV-version only)
- Base plate included, angle bracket included (5SV-version only)

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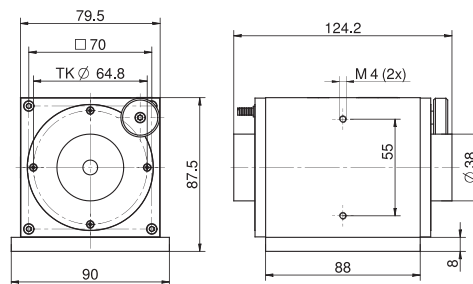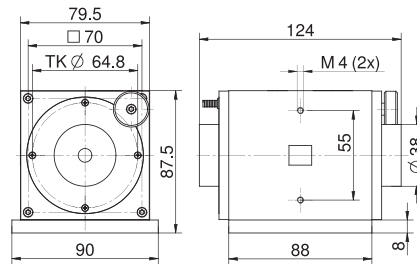
Tunable Isolator with 8 mm Aperture, SI Series

FI-x/y-8SI

- Continuous adjustment for wavelength without movement of optical components
- Tunable with maximum transmission and isolation over the complete wavelength range
- Isolation better than 30 dB, typically 38-42 dB over the entire wavelength range
- Custom isolation values on request
- TGG crystal
- Access to blocked beam
- Mounting: via two M4 threaded holes at the bottom side and the back side, 55 mm separation, or via base plate
- Base plate (88x90x8 mm³) included

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A precise mechanics enables a continuous wavelength adjustment. Without movement of the optics a broad wavelength range is realized.

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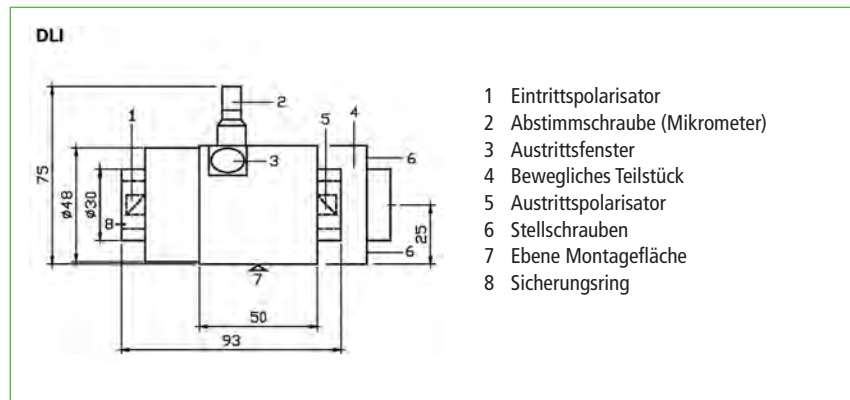
Two Stage Faraday Isolators, Technical Overview

FI-x-5TI and FI-x-5TV

Diode lasers are extremely sensitive to reflected radiation. Standard Faraday isolators typically achieve between 30dB and 40 dB isolation, which in some cases is not sufficient to suppress undesirable feedback.

Our two stage Faraday isolators were developed for the special requirements of diode lasers and square the standard isolation of single stage Faraday isolators. At the heart of this development is the use of two coupled isolator stages together with the best polarizers available on the market.

This configuration combines the exit polarizer of the first stage with the entry polarizer of the second stage to form one central polarizer.

Arranging the polarity of the two magnets to be opposite to each other results in two benefits: The polarization direction of the transmitted light remains unchanged in the transmission direction and the effect of both magnetic fields is enhanced. Therefore this configuration also leads to a more compact isolator and a reduction of the optical path length which in turn enhances the optical quality of the Faraday isolator.

All optical surfaces are antireflection coated and the surfaces normal to the beam axis are tilted. The polarizers are mounted in a way that allows easy cleaning of the external optical surfaces. This guarantees that the isolation is not reduced by residual reflections and scattering from the isolator.

Based on this special design a guaranteed 60 dB isolation at the design wavelength, and within the tuning range makes our two stage Faraday isolators the best on the market.

A precise wavelength adjustment is possible by rotating the central polarizer with a micrometer set screw. The blocked radiation is deflected out of the isolator at 90° with respect to the beam axis. It is not absorbed by the interior of the isolator, but is available at the side surfaces of the polarizer and the exit window.

DLI Injection Locking

The DLI injection version revolves this operating mode and uses the exit window for in-coupling of the seed laser for injection locking while decoupling efficiently the master and the slave laser from each other at the same time. Like this stable mode locking (e.g. of Ti: Sapphire lasers) is simplified.

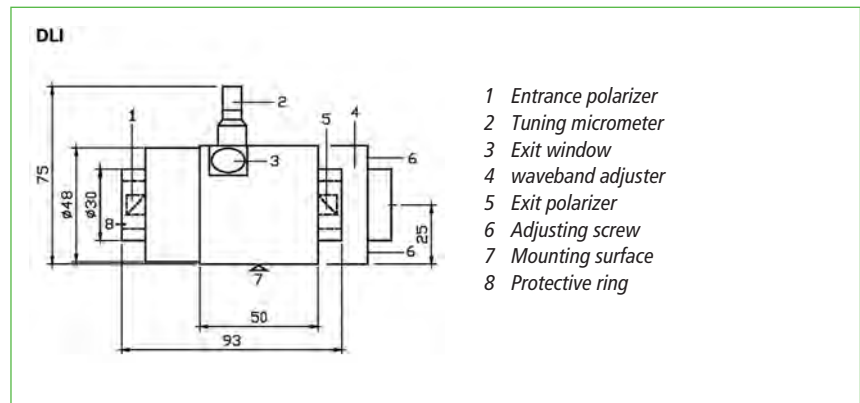
Applications

All two stage isolators are typically used to improve the power and frequency stability of diode lasers used in spectroscopy, interferometry and precision control as well as in alignment applications. Since the output polarization and the beam position are conserved for all two stage Faraday isolators, the influence of the smallest feedback effects on the laser can be quantitatively examined.

DLI, Overview

The isolators of the DLI-series were developed for the special requirements of diode lasers in the visible spectrum and combine the outstanding isolation of a two stage isolator with the flexibility of a tunable isolator.

The DLI isolators are easily integrated into an existing setup and can be adjusted to match any wavelength without changing the laser polarization or displacing the laser beam. The isolators can be coarsely tuned by altering the effective magnetic field in the two isolator stages.

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
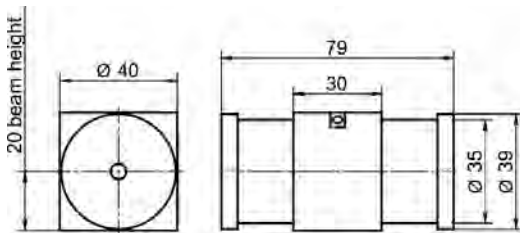
Two-Stage Faraday Isolators, tunable, TVC/TIC Series

FI-x-TV / FI-x-TI

- Tunable with maximum isolation over the complete wavelength range
- Two coupled isolator stages in series
- Especially high isolation > 60 dB
- TGG crystal
- Customized central wavelength on request
- Mounting TVC-Version: via two M4 threaded holes on the bottom side and at the back side, 30 mm separation
- Mounting TIC-Version: via 8xM2 threaded holes on the bottom side and on the back side, 22.5 mm separation

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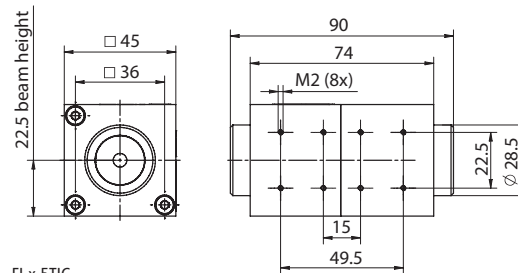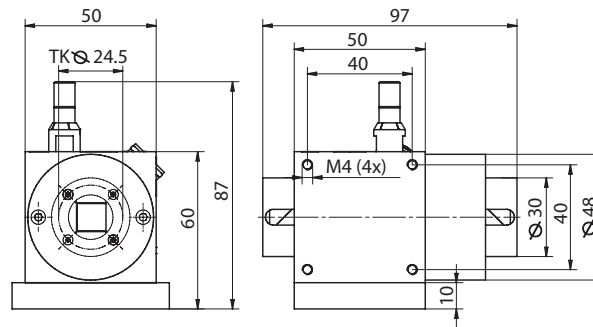
Two-Stage Faraday Isolators, tunable, DLI Series

Tunable Diode Laser Isolators DLI

- Tunable with maximum isolation and maximum transmittance over the complete wavelength range
- Two coupled isolator stages in series
- Especially high isolation ≥ 60 dB
- TGG crystal
- Input polarization = output polarization
- Individually calibrated adjustment curve supplied with each isolator
- Aperture $\varnothing 5$ mm
- Mounting: via four M4 threaded holes at the bottom side and at the back side, 40x40 mm separation, or via base plate
- Base plate (50x60x10 mm³) included
- Access to blocked beam via window

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An easy integration of DLI isolators is possible. They can be adjusted without changing laser polarization or beam position. The outer window can be used for injection locking.

Zweistufige Isolatoren, durchstimbar, DLI-Reihe Two-Stage Faraday Isolators, tunable, DLI Series

Produktbezeichnung	Isolation, garantiert (dB)	Transmission bei den Designwellenlängen (%)	Einstellbereich (typisch) (nm)	Abmessungen (mm)	Zerstörschwelle @ 1.2 ps Pulse (J/cm ²)	Zerstörschwelle @ 8 ns Pulse (J/cm ²)	Part No.
Item Title	Isolation, guaranteed (dB)	Transmission at design wavelength (%)	Tuning range typical (nm)	Dimensions (mm)	Damage threshold @ 1.2 ps pulses (J/cm ²)	Damage threshold @ 8 ns pulses (J/cm ²)	Part No.
DLI 1	≥ 60	≥ 80	745 - 890	50x50x97	0.2 @ 890 nm	0.9 @ 890 nm	84501003000
DLI 2	≥ 60	≥ 80	610 - 700	50x50x97	0.1 @ 700 nm	0.7 @ 700 nm	84501002000
DLI 3	≥ 60	≥ 80	650 - 760	50x50x97	0.2 @ 760 nm	0.8 @ 760 nm	84501001000

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Our quality criteria:

- *Best possible extinction ratio for each crystal*
- *High transmission*
- *Patented isolation system minimizes piezoelectric oscillation for exceptionally precise switching operations (optional)*

Laser Modulators & Pockels Cells

In the following section electro-optical modulators are divided into modulators (for applications outside of laser cavity) and Pockels cells (for applications within laser cavity).

You can choose from a large selection of crystals for a variety of applications, apertures and laser powers, covering the entire wavelength range from 250 nm to 3 μm . The consistently high Qioptiq quality and incomparable value of our products is assured by a combination of our many years of experience, an intelligent design, modern engineering with computer simulations and sophisticated processing.

In addition we offer a broad range of fast high-voltage drivers.

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Special Features:

If technically feasible, we can modify any product for wavelengths in the 250 nm to 3 μm range on request..

Ideal Areas:

Phase and intensity modulation; Q-switching; pulse picking.

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Pockels Cells

<i>Technical Information</i>	544
<i>Product Overview</i>	546
<i>KD*P Pockels Cells LM Series</i>	547
<i>KD*P Pockels Cells CPC Series</i>	548
<i>KD*P Pockels Cells CIQS Series</i>	549
<i>KD*P Pockels Cells SPC 4 Series</i>	550
<i>KD*P Double Pockels Cells DPZ Series</i>	551
<i>KD*P Brewster Pockels Cell BPC 8</i>	552
<i>BBO Pockels Cells BBPC Series</i>	552
<i>BBO Double Pockels Cells DBBPC Series</i>	553
<i>BBO Pockels Cells for High Switching Rates DBBPC HR Series</i>	554
<i>RTPC Pockels Cells Series</i>	555
<i>Pockels Cells Positioner</i>	556

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Introduction and Technical Overview

Electro-optical crystals are characterized by their ability to change optical path length in function of an applied external voltage. This change depends on the direction of polarization of the irradiated light. At $N/2$ voltage, the path length difference of orthogonally polarized beams is just half of the wavelength. With a suitable orientation of the crystals, the polarization direction of the irradiated light is rotated 90° : in this state the light is extinguished by a polarizer. Varying the applied voltage allows quick modulation of the laser beam intensity. The performance of an electro-optic modulator can be understood very simply as that of a retardation plate with electrically adjustable retardation.

Series LM 0202 Modulators use the transverse electro-optical effect: the direction of the light beam and electric field are orthogonal. In this configuration, long crystals with a small cross section have a low half wave voltage

Since most of the electro-optical crystals operate with a strong background of natural birefringence, a compensation scheme is used. Each modulator in the LM 0202 series has four crystals as a matched ensemble. These crystals are fabricated with deviations in length less than 100 nm. The crystals are operated optically in series and electrically parallel.

The crystal orientation of the LM 0202 and LM 0202P modulators has been optimized to minimize the retardation caused by natural birefringence. Just as in an ordinary retardation plate, the polarization of the laser beam has to be adjusted at 45° to the optical axis in order to achieve a proper 90° rotation.

If the laser beam is polarized in the direction of the optical axis, no polarization rotation, but pure phase retardation will occur. In principle this allows the user to operate the LM 0202 modulator as a phase modulator. In this configuration, optimized for minimum background retardation, two of the four crystals are electro-optically active for phase modulation. A special model, LM 0202 PHAS, is available with a crystal configuration that uses all four crystals for phase modulation.

The PM 25 Phase Modulator, is a Brewster modulator of high optical quality and should be used for loss sensitive applications, especially intracavity modulation. Mounting the modulator in the resonator is simple, as there is no beam deviation or displacement.

For all modulators, the beam deflection and beam displacement caused by the natural birefringence is eliminated by using the crystals in compensation configuration.

Electro-optic modulators generally require linearly polarized laser light. If the laser light is not sufficiently polarized by itself, an additional polarizer must be used.

The Intensity Modulator LM 0202 P has an integrated polarizer that is used as an analyzer.

The modulator voltage input plugs are isolated from the housing and directly connected to the crystals. A change of the laser intensity can be observed when the applied voltage is changed. By subsequently adjusting voltage and rotation, an extinction better than 250:1 can be achieved. Selected models with better extinction ratios are available on request.

Operating an electro-optical modulator between crossed, or parallel, polarizers yields an intensity variation given by the following formula:

$$I = I_0 \cdot \sin^2 (U/U_{N/2} \cdot \pi/2)$$

$U_{N/2}$ - half wave voltage
 I_0 - input intensity
 U - signal voltage

It has been assumed that the appropriate offset voltage has been applied for maximum extinction. The offset voltage causes a shift of the intensity curve over the voltage. The half wave voltage is proportionally to the wavelength λ , to the crystal thickness d and in reverse proportional to the crystal length l :

$$U_{N/2} = \frac{\lambda \cdot d}{n_0^3 r_{63} l}$$

Here n_0 is the refractive index of the ordinary beam and r_{63} the electro-optical coefficient of the crystal.

In many cases it is advantageous to select an offset voltage such that the first order intensity varies linearly with voltage. This is achieved by setting the offset voltage to the value required for maximum extinction minus $\frac{1}{2} \cdot U_{N/2}$

The LM 0202 series modulators are hermetically sealed. They can be operated at pressures from 100 mbar to 1500 mbar and at a temperature range between 0°C to 50°C .

Standard models are designed for horizontal operation. Modulators for vertical use are available by request. The modulator windows are easily cleaned with a mild organic solvent.

Applications

The series LM 0202 or LM 13 electro-optical modulators are typically used when intensity, power, phase or polarization state modulation is required. The devices are ideal for continuous or pulsed laser applications. Standard models, in many configurations, are available for wavelength ranges or for definite wavelengths between 250 to 1100 nm and operation up to 3000 nm is possible with special crystals.

The modulators are typically used with diode lasers, solid state lasers, ion lasers, gas lasers or white light lasers.

These devices are being used in the fields of laser scanning microscopes, stereo lithography, laser projection, optical storage, printing, research and development and communication engineering in the laser industry.

The PM 25 and PM-C-BB series are typically used for fast intra-cavity phase modulation. Therefore very fast control loops, with high feedback gain for frequency and phase stabilization, can be constructed for precision lasers.

Selection Criteria

The required wavelength and aperture are determined based on the existing laser system. Very high laser power, in the multiwatt range, requires a large aperture. Laser lines in the short wave spectral region can work easily with modulators having low electro-optical sensitivity: this gives rise to advantages in bandwidth and size. A Brewster modulator of high optical quality should be used for loss sensitive applications, especially intracavity modulation.

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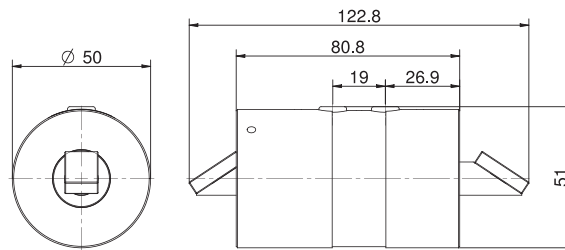
Phase Modulator PM 25

- Two crystals at Brewster angle in order of compensation
- With Brewster windows
- Very high transmission
- Connectors: 4 mm banana plugs
- Different versions for wavelength ranges between 300 and 1100 nm
- Wavefront Distortion: $< \lambda/10$ at 633 nm
- Bandwidth (3 dB): 100 MHz
- Capacitance: 30 pF
- Max. continuous voltage: 1500 V
- Operating temperature: 10-45 °C
- Weight: 500 g (17.6 oz) approx.
- Please specify the wavelength or wavelength range and laser parameters when ordering
- Modifications and custom specifications on request

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
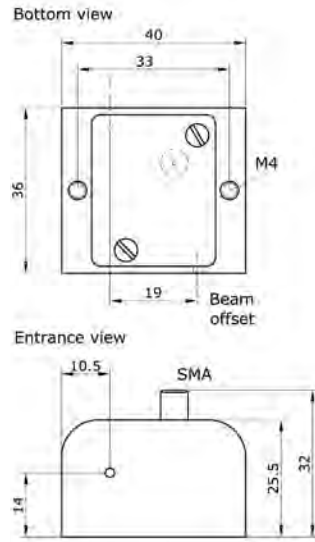
Phase Modulator PM-C-BB

- Brewster-cut MgO-LiNbO₃ crystal
- Broad wavelength range 450-3000 nm (except LiNbO₃ absorption at 2.82-2.84 μm)
- High photorefractive damage threshold > 100 W/mm² @1064 nm (cw operation, wavelength dependent)
- High transmission
- Small residual amplitude modulation
- Compact design
- Capacitance: 10 pF
- PM-C-BB (T) with built-in active temperature stabilization (< 10 mK)
- Wavefront distortion: < λ/4 at 633 nm
- Bandwidth: DC-500 MHz (> 10 MHz resonance-free)
- Connector: 1 x SMA

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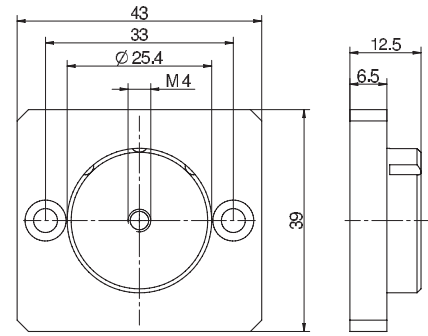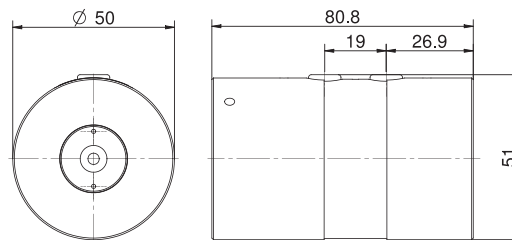
Laser Modulators LM 13

- Different versions: universal modulator, intensity modulator (P) with thin film polarizer, phase modulator (PHAS)
- With 2 crystals in order of compensation
- Connectors: 4 mm banana plugs
- Different versions for wavelength ranges between 250 and 1100 nm
- Extinction: > 250:1 (VIS, IR) or > 100:1 (UV) (measured at continuous wave between crossed polarizers)
- Wavefront Distortion $< \lambda/4$ at 633 nm
- Bandwidth (3 dB): 100 MHz
- Capacitance: 46 pF
- Max. continuous voltage: 800 V
- Operating temperature: 10-45 °C
- Weight: 500 g (28.2 oz) approx.
- Please specify the wavelength or wavelength range and laser parameters when ordering
- Modulators series LM 13 are also available with LiTaO₃ crystals for a broader wavelength range
- Modifications and custom specifications on request

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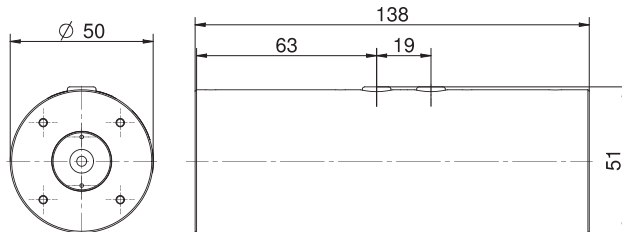
Laser Modulators LM 0202

- Different versions: universal modulator, intensity modulator (P) with thin film polarizer, phase modulator (PHAS)
- With 4 crystals in order of compensation
- Connectors: 4 mm banana plugs
- Different versions for wavelength ranges between 250 and 1100 nm
- Extinction: > 250:1 (VIS, IR) or > 100:1 (UV) (measured at continuous wave between crossed polarizers)
- Wavefront Distortion $< \lambda/4$ at 633 nm
- Bandwidth (3 dB): 100 MHz
- Capacitance: 82 pF
- Max. continuous voltage: 800 V
- Operating temperature: 10-45 °C
- Weight: 800 g (17.6 oz) approx.
- Please specify the wavelength or wavelength range and laser parameters when ordering
- Modulators series LM 0202 are also available with LiTaO₃ crystals for a broader wavelength range
- Modulators series LM 0202 are also available with ADP crystals for an even more precise performance
- Modifications and custom specifications on request

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Digital Pulse Amplifier DIV 20

- Digital pulse amplifier for all LM0202 and LM13 laser modulators with half wave voltage up to 590 V
- Generates rectangular voltage pulses of variable pulse length and adjustable amplitude
- Control via 5 V TTL signal

Specifications:

Signal voltage ¹⁾²⁾: 30-590 V
Rise/fall time (10-90 %) ³⁾: < 28 ns, (typ. 20-25 ns at 350 V, typ. 23-27 ns at 590 V)

Repetition rate ⁴⁾: 1 to 20 MHz
Offset-voltage ^{1) 2)}: 0-590 V
External control via Sub-D25p-female-port
Line voltage: 90-264 V
Line frequency: 47-63 Hz

¹⁾ Relative to ground

²⁾ This voltage can be set manually or externally by a control voltage from 0 to +10 V (input impedance 5 kΩ) with a speed of <100 ms from 70 V to 480 V (rise) or from 480 V to 70 V (fall) respectively

³⁾ The optical rise/fall time is achieved with a modulator LM 0202, connected via the standard connection cable (L = 80 cm)

⁴⁾ Typical maximum signal voltage at repetition rate 5 MHz 300 V, at 3 MHz 400 V and at 1 MHz 590 V

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SINE-Amplifier for Phase Modulators

- Compact design
- Can be used with PM-C-BB, PM 25, LM 13 and LM 0202
- Large modulation bandwidth
- High output voltage
- Cost effective
- Modulator cable and adapters included
- Power-supply included (with universal power supply and snap-in adapters for EU, US, GB, AUS)
- Input waveform: SINE-Wave
- Input voltage: max. +13 dBm
- Bandwidth with LM 0202: appr. 4-7 MHz
- Bandwidth with PM-C-BB: appr. 5-12 MHz
- Max. output voltage with LM 0202: 200 V_{pp} @ 5-7 MHz frequency
- Max. output voltage with PM-C-BB appr.: > 500 V_{pp} @ 8-12 MHz frequency
- Supply-voltage (power-supply incl.): +12 V DC
- Dimensions: 115x65x70 mm³

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Technical Information

The Electro-Optic Effect

The linear Electro-Optic effect, also known as the Pockels effect, describes the variation of the refractive index of an optical medium under the influence of an external electrical field. In this case certain crystals become birefringent in the direction of the optical axis which is isotropic without an applied voltage.

When linearly polarized light propagates along the direction of the optical axis of the crystal, its state of polarization remains unchanged as long as no voltage is applied. When a voltage is applied, the light exits the crystal in a state of polarization which is in general elliptical.

This way phase plates can be realized in analogy to conventional polarization optics. Phase plates introduce a phase shift between the ordinary and the extraordinary beam. Unlike conventional optics, the magnitude of the phase shift can be adjusted with an externally applied voltage and a $\lambda/4$ or $\lambda/2$ retardation can be achieved at a given wavelength. This presupposes that the plane of polarization of the incident light bisects the right angle between the axes which have been electrically induced. In the longitudinal Pockels effect the direction of the light beam is parallel to the direction of the electric field. In the transverse Pockels cell they are perpendicular to each other. The most common application of the Pockels cell is the switching of the quality factor of a laser cavity.

Q-Switching

Laser activity begins when the threshold condition is met: the optical amplification for one round trip in the laser resonator is greater than the losses (output coupling, diffraction, absorption, scattering). The laser continues emitting until either the stored energy is exhausted, or the input from the pump source stops. Only a fraction of the storage capacity is effectively used in the operating mode. If it were possible to block the laser action long enough to store a maximum energy, then this energy could be released in a very short time period to form a giant light pulse.

A method to accomplish this is called Q-switching. The resonator quality, which represents the losses in the resonator, is kept low until the maximum energy is stored. A rapid increase of the resonator quality then takes the laser high above threshold, and the

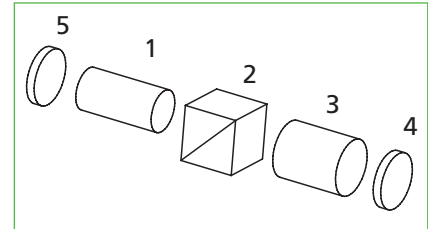
stored energy can be released in a very short time. The resonator quality can be controlled as a function of time in a number of ways. In particular, deep modulation of the resonator quality is possible with components that influence the state of polarization of the light. Rotating the polarization plane of linearly polarized light by 90° , the light can be guided out of the laser cavity by a polarizer.

The modulation depth, apart from the homogeneity of the 90° rotation, is only determined by the extinction ratio of the polarizer. The linear electro-optical (Pockels) effect plays a predominant role compared to the quadratic electro-optical (Kerr) effect. Typical electro-optic Q-switches operate in a so called $\lambda/4$ mode.

a) Off-Q-Switching


Light emitted by the laser rod (1) is linearly polarized by the polarizer (2). If a $\lambda/4$ voltage is applied to the Pockels cell (3), then on exit, the light is circularly polarized. After reflection from the resonator mirror (4) and a further passage through the Pockels cell, the light is linearly polarized again, but the plane of polarization has been rotated by 90° . The light is deflected out of the resonator at the polarizer. Hence the resonator quality is low and the laser does not start to oscillate. At the moment the maximum storage capacity of the active medium has been reached, the voltage of the Pockels cell is turned off very rapidly; the resonator quality increases immediately and a very short laser pulse is emitted. The use of a polarizer can be omitted for active materials which show polarization dependent amplification (eg. Nd:YAlO₃, Alexandrite, Ruby, etc.).

Off-Q-Switching

b) On-Q-Switching

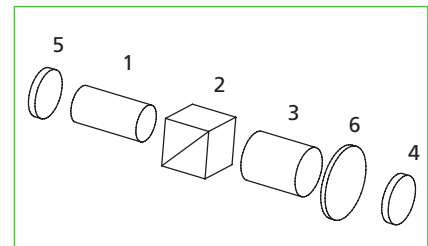Unlike off Q-switching, a $\lambda/4$ plate (6) is used between the Pockels cell (3) and the resonator mirror (4). If no voltage is applied to the Pockels cell the laser resonator is blocked: no laser action takes place. A voltage pulse then compensates for the waveplate, thus opens the resonator and permits the emission of laser light.

On-Q-Switching

Pulse Picking

Typically Femto-Second-Lasers emit pulses with a repetition rate of several 10 MHz. However many applications like regenerative amplifying require slower repetition rates. Here a Pockels cell can be used as an optical switch: by applying ultra fast and precisely timed $\lambda/2$ -voltage pulses on the Pockels cell, the polarization of the Laser light can be controlled pulse wise. Thus, combined with a polarizer the Pockels cell works as an optical gate.

Selection Criteria

The selection of the correct Q-switch for a given application is determined by the excitation of the laser, the required pulse parameters, the switching voltage, the switching speed of the Pockels cell, the wavelength and the polarization state of the light.

Type of Excitation

Basically, both off and on Q-switching are equivalent in physical terms for both cw and for pulse pumped lasers. On-Q-switching is, however, recommended in cw operation because a high voltage pulse and not a rapid high voltage switch-off is necessary to generate a laser pulse. This method also extends the life time of the cell. Over a long period of time, the continuous application of a high voltage, as well as $\lambda/2$ -voltage applications would lead to electrochemical degradation effects in the KD*P crystal. We advise the use of an On-Q-switching driver.

Off Q-switching is more advantageous for lasers stimulated with flash lamps because the $\lambda/4$ plate is not required. In order to prevent the electrochemical degradation of the KD*P crystal in the off Q-switching mode we recommend a trigger scheme in which the high voltage is turned off between the flashlamp pulses and turned on to close the laser cavity before the onset of the pump pulse.

The cell CPC and SPC series are recommended for diode pumped solid state lasers. These cells are ultra compact and will operate in a short length resonator: this is necessary to achieve very short laser pulses.

Pulse Parameters

The series LM n IM, and LM n SG cells are recommended for lasers with a power density of up to 2 GW/cm². The LM n and LM n SG cells are used for lasers with very high amplification. The SG cells with sol-gel technology have the same transmission as the immersion cells and both are typically used when a higher transmission is required.

Brewster Pockels cells are recommended for lasers with low amplification, such as Alexandrite lasers. The passive resonator losses are minimal due to a high transmission of 99 %.

The CPC and SPC series cells are suitable for small, compact lasers and especially for OEM applications. They are available as Sol-Gel cells and immersion cells.

The level of Deuterium content in a KD*P crystal influences the spectral position of the infrared edge. The higher the deuterium level the further the absorption edge is shifted into the infrared spectral region: for Nd: YAG at 1064 nm, the laser absorption decreases. Crystals, which are deuterated to > 98 %, are available for lasers with a high repetition rate or a high average output power.

Pockels Cell Switching Voltage

Using double Pockels cells can reduce the switching voltage by 50%. This is achieved by switching two crystals electrically in parallel and optically in series. The damage threshold is very high and the cells are mainly used outside the resonator.

Electro optic material

The selection of the electro-optic material depends on its transmission range. Further the laser parameters and the application as well have to be taken into account.

For wavelengths from 0.25 μm to 1.1 μm , longitudinal Pockels cells made of KD*P and a deuterium content of 95 % should be considered. If the deuterium content is higher the absorption edge of the material is shifted further into the infrared. KD*P crystal cells with a deuterium content > 98 % can be used up to 1.3 μm .

KD*P can be grown with high optical uniformity and is therefore recommended for large apertures.

The spectral window of BBO also ranges from 0.25 μm to 1.3 μm . In addition BBO crystals provides a low dielectric constant and a low absorption. Therefore BBO is recommended for Lasers with high repetition rate and high average powers.

RTP, with an optical bandwidth from 0.5 μm up to 1.5 μm , combines low switching voltage and high laser induced damage threshold. Together with its relative insensitivity for Piezo effects RTP is best suited for precise switching in high repetition rate lasers with super fast voltage drivers.

Suppression of Piezo effects

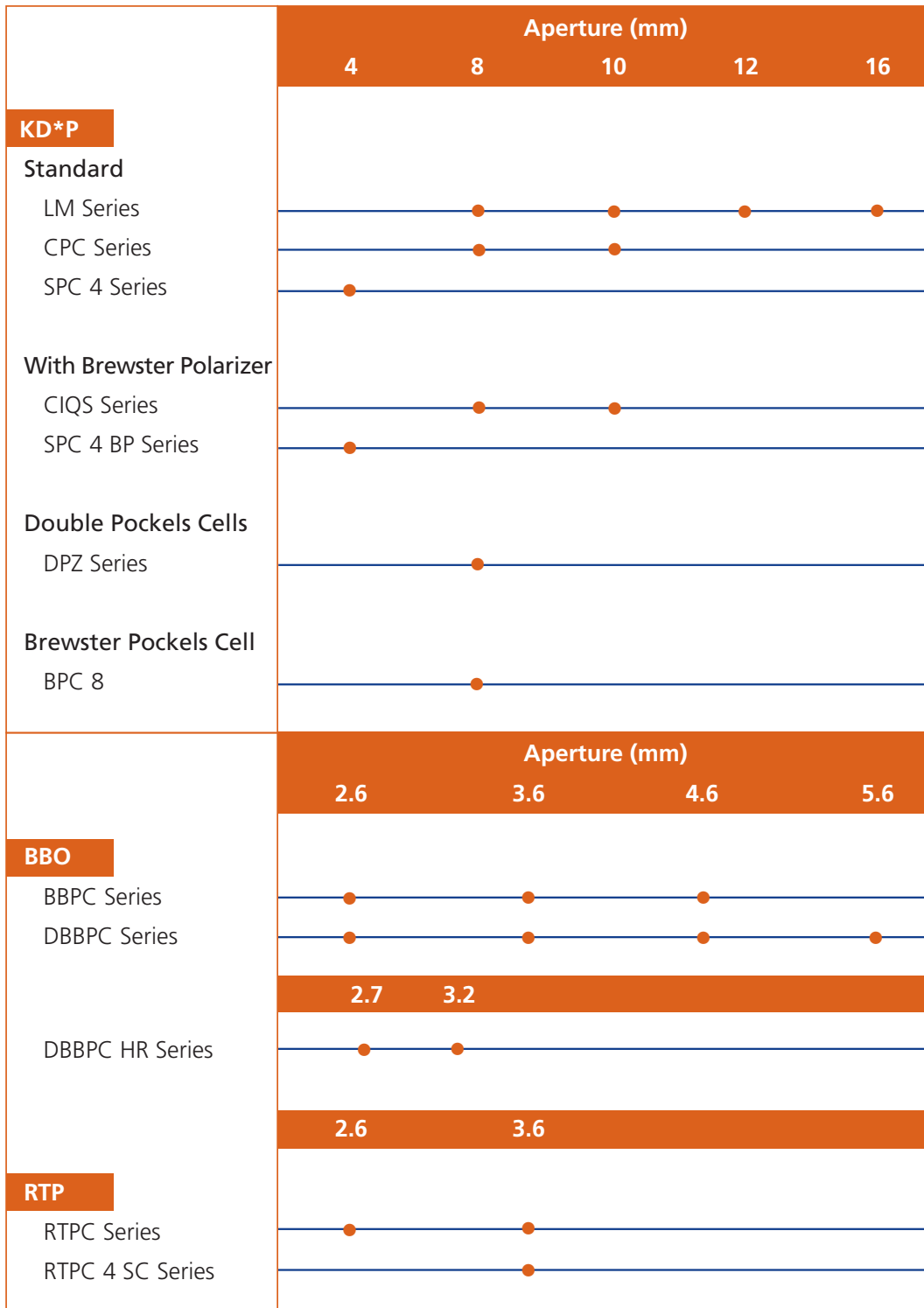
Like any other insulating material electro optical crystals show Piezo effects when high voltage is applied. The extend of the Piezo ringing depends on the electro optic material and usually its effect on the extinction ratio is negligible when used for Q-switching. However for pulse picking applications, which require highly precise switching behaviour, LINOS offers specially Piezo damped Pockels cells which suppress these ringing effects efficiently.

State of Polarization

The CIQS series cells are supplied with an integrated polarizer and an optical waveplate: The alignment of the Pockels cell relative to the polarizer thus can be omitted. The rotational position of the cell relative to the resonator axis can be chosen at will. However, should the polarization state of the light in the resonator be determined by other components, such as anisotropic amplification of the laser crystal or Brewster surfaces of the laser rod, then the rotational position of the cell will be determined by these factors. Thin film polarizers are used and the substrate is mounted at the Brewster angle. A parallel beam displacement of 1 mm results from this configuration and can be compensated by adjusting the resonator easily.

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

KD*P Pockels Cells LM Series

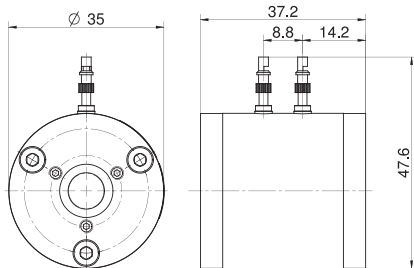
- KD*P-based pockels cells
- High crystal deuteration (typical): > 98%
- Wavefront Distortion: < $\lambda/4$
- Damage threshold: > 2 GW/cm² at 1064 nm, 12 ns, 1 Hz (typical)
- Available as immersion (IM) or SolGel (SG) version
- All order numbers valid for 1064 nm
- Other specifications on request
- Please state the applied wavelength when ordering.

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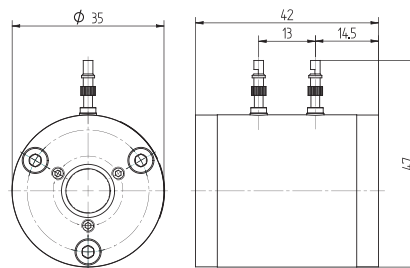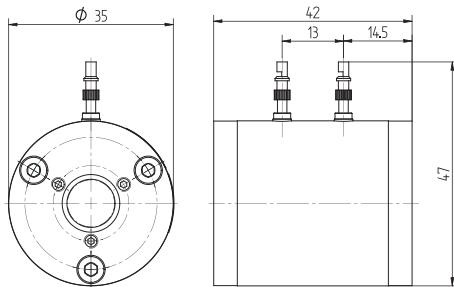
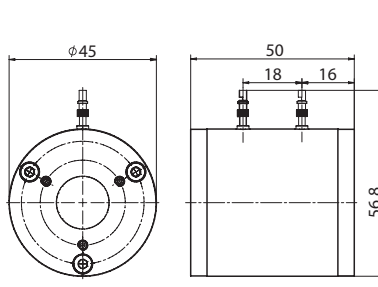
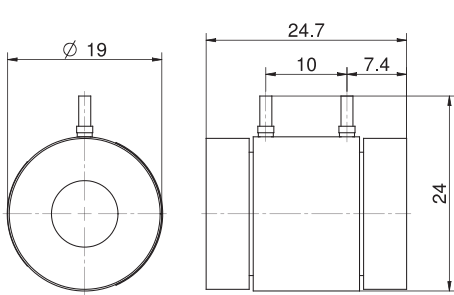
KD*P Pockels Cells CPC Series

- KD*P-based pockels cells
- Compact design for OEM applications
- High crystal deuteration (typical): > 98%
- Wavefront Distortion: < $\lambda/4$
- Damage threshold: 2 GW/cm² at 1064 nm, 12 ns, 1 Hz (typical, not guaranteed)
- Available as immersion (IM) or SolGel (SG) version
- Other specifications on request
- Please state the applied wavelength when ordering.

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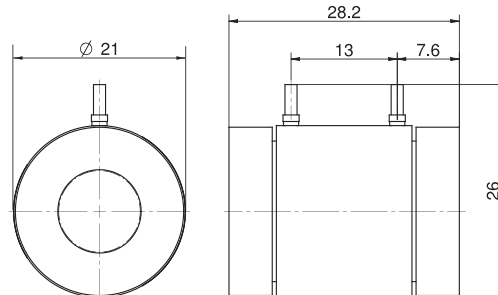
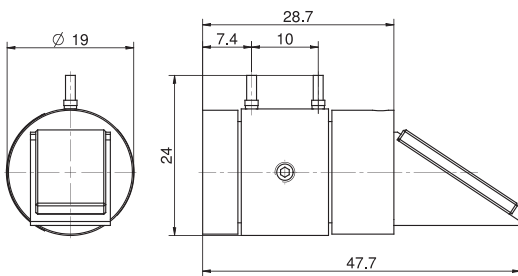
KD*P Pockels Cells CIQS Series


- KD*P-based pockels cells
 - With integrated, pre-adjusted Brewster polarizer
 - Compact design for OEM applications
 - High crystal deuteration (typical): > 98%
 - Wavefront Distortion: < $\lambda/4$
 - Damage threshold: 1.5 GW/cm² at 1064 nm, 12 ns, 1 Hz (typical, not guaranteed)
- Available as immersion (IM) or SolGel (SG) version
 - All order numbers valid for 1064 nm
 - Other specifications upon request
 - Please state the applied wavelength when ordering.
 - Versions with integrated $\lambda/4$ -waveplate available

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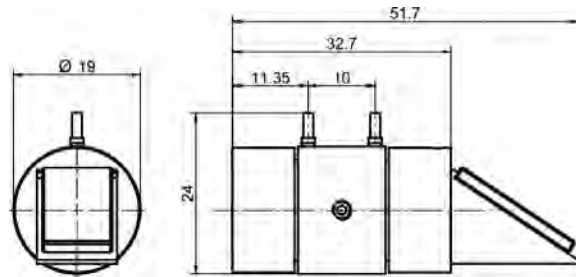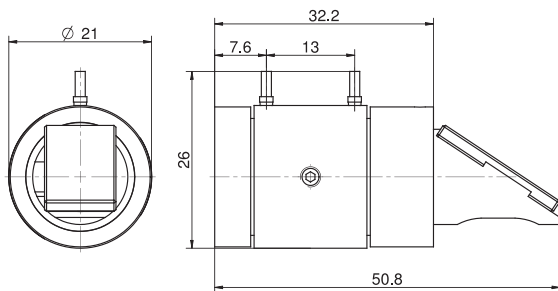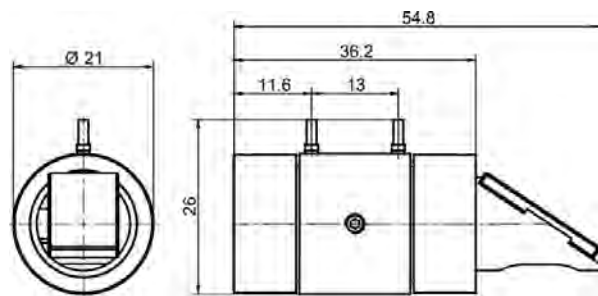
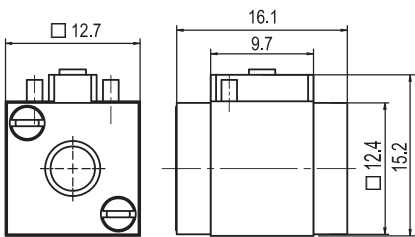
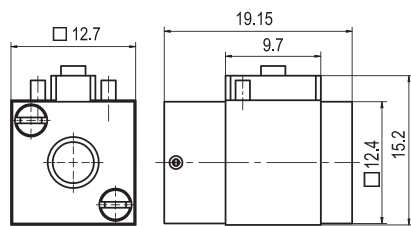
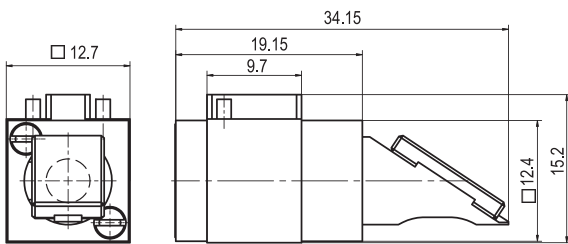
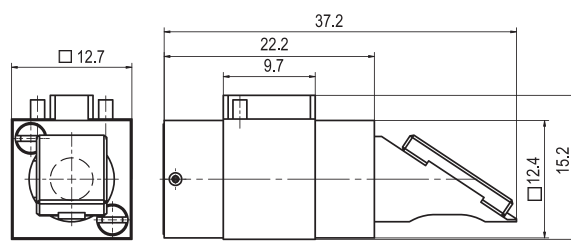
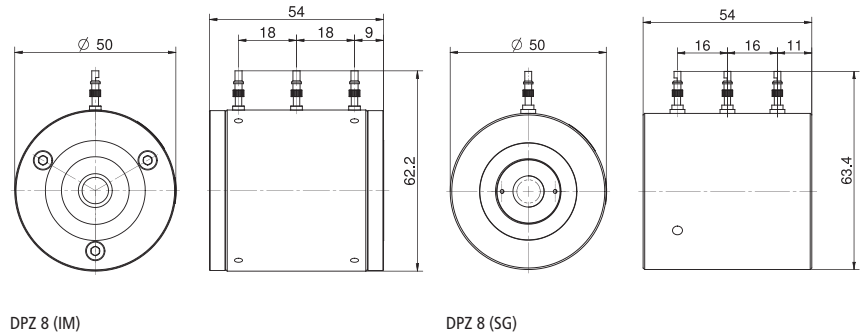
KD*P Double Pockels Cells DPZ Series

- KD*P-based pockels cells
- Two crystals in series reduce the switching voltage by 50%
- High crystal deuteration (typical): > 98%
- For shutter and $\lambda/2$ -applications
- Damage threshold: 2 GW/cm² at 1064 nm, 12 ns, 1 Hz (typical)
- Available as immersion (IM) or SolGel (SG) version
- $\lambda/4$ -voltage: 1.6 kV at 1064 nm, 20°C
- All order numbers valid for 1064 nm
- Other specifications on request
- Please state the applied wavelength when ordering.


KD*P Doppel-Pockelszelle der DPZ-Serie - 1064 nm KD*P Double Pockels Cells DPZ Series - 1064 nm

Produktbezeichnung	Apertur Ø (mm)	Transmission, typ. (%)	Löschungsverhältnis (spannungsfrei)	$\lambda/2$ -Spannung at 1064 nm, 20°C (kV)	Kapazität (pF)	Part No.
Item Title	Aperture Ø (mm)	Transmittance, typ. (%)	Extinction ratio (voltage-free)	$\lambda/2$ -voltage at 1064 nm, 20°C (kV)	Capacitance (pF)	Part No.
DPZ 8 IM	8	95	> 1000:1	3.2 +15% / - 10%	8	84503042000
DPZ 8 SG	7.5	95	> 1000:1	3.2 +15% / - 10%	8	84503043005

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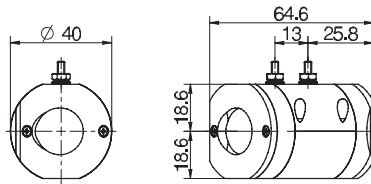
KD*P Brewster Pockels Cell BPC 8

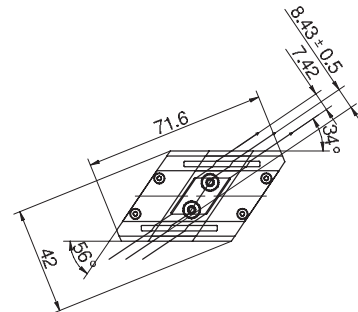- KD*P-based pockels cells
- High crystal deuteration (typical): > 98%
- Crystal with Brewster angle cut
- High transmittance for lasers with low amplification
- Mounting via ENSAT-SBI M4 thread insert at the bottom side
- Beam offset: 8.4 mm
- Wavefront Distortion: < $\lambda/4$
- Damage threshold: > 500 MW/cm² at 1064 nm, 10 ns, 1 Hz (typical)
- Other specifications on request.
- Please state the applied wavelength when ordering.

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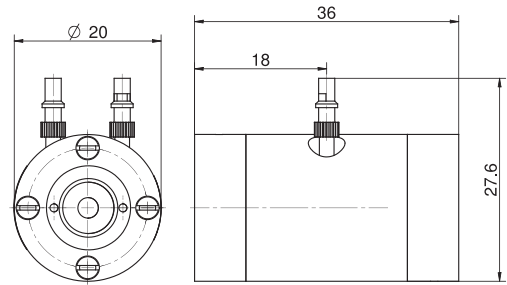
BBO Pockels Cells BBPC Series

- BBO-based pockels cells
- Suited for Q-switch applications with high repetition rates
- Wavefront Distortion: < $\lambda/4$
- Damage threshold: > 4 J/cm² @ 74 ps, 1064 nm
- Optionally available with integrated Brewster polarizer: BBPC n BP
- Optionally available with integrated $\lambda/4$ waveplate: BBPC n WP
- Optionally available with piezo attenuator: BBPC n pp
- All order numbers valid for 1064 nm
- Other specifications on request.
- Please state the applied wavelength when ordering.

BBPC

BBO-Pockelszellen der BBPC-Serie - 1064 nm
BBO Pockels Cells BBPC Series - 1064 nm

Produkt-bezeichnung	Apertur Ø (mm)	Transmission, typ. (%)	Löschungsverhältnis (spannungsfrei)	λ/4-Spannung DC at 1064 nm, 20°C (kV)	Kapazität (pF)	Part No.
<i>Item Title</i>	<i>Aperture Ø (mm)</i>	<i>Transmittance, typ. (%)</i>	<i>Extinction ratio (voltage-free)</i>	<i>λ/4-voltage DC at 1064 nm, 20°C (kV)</i>	<i>Capacitance (pF)</i>	<i>Part No.</i>
BBPC 3	2.6	98	>1000:1	3.6 ± 15%	4	84503083012
BBPC 4	3.6	98	>1000:1	4.8 ± 15%	4	84503083008
BBPC 5	4.6	98	>1000:1	6.0 ± 15%	4	84503083020

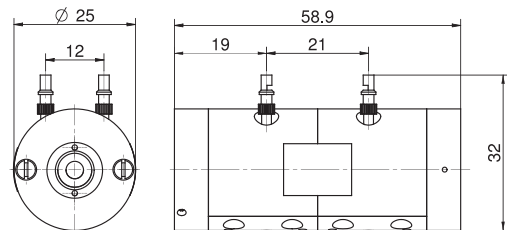
BBO-Doppel-Pockelszellen der DBBPC-Serie

BBO Double Pockels Cells DBBPC Series

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- BBO-based double pockels cells
- Two crystals in series reduce the switching voltage by 50%
- With piezo attenuator
- Suited for Q-switch applications with high repetition rates, pulse picking and regenerative amplifiers
- Damage threshold: > 4 J/cm² @ 74 ps, 1064 nm
- All order numbers valid for 1064 nm
- Other specifications on request
- Please state the applied wavelength when ordering.

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
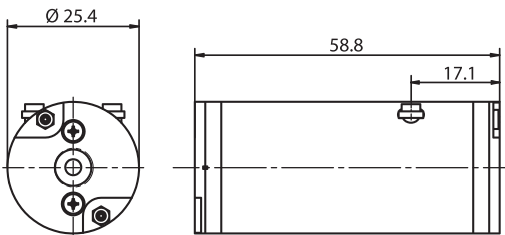
BBO Pockels Cells for High Switching Rates DBBPC HR Series

The new LINOS Double BBO High-Repetition Pockels Cells DBBPC HR from Qioptiq enable the fastest possible switching rates available today with > 1.3 MHz performance. Together with the low absorbing BBO crystals the DBBPC HR Pockels cells are perfectly suited for the next generation of fast high power regenerative amplifiers and pulse pickers.

- BBO Pockels Cell with 2 crystals in optical series
 - For high switching rate applications up to > 1.3 MHz
 - For high average power applications
 - Design wavelength 1030 nm (other wavelengths on request)
 - Damage threshold @ 1064 nm, 74 ps: > 4 J/cm²
 - Reduced switching voltage
 - Suitable for quarter- and half-wave mode applications
 - For Q-Switching, cavity dumping, regenerative amplifiers and pulse pickers
- All order numbers valid for 1030 nm
 - Other specifications on request
 - Please state the applied wavelength when ordering.
 - This Product is export controlled according to regulation (EC) No. 428/2009, list item 6A005. Shipping is subject to the existence of a valid export license, if necessary.

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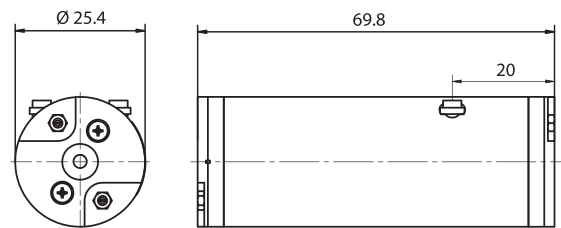
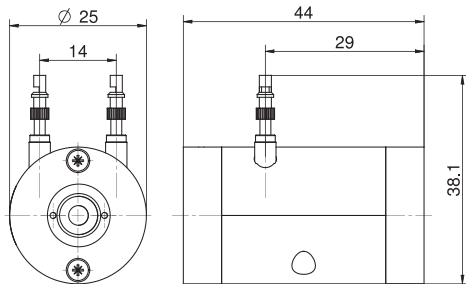
RTPC Pockels Cells Series

- RTP-based pockels cells
- Suited for Q-switch applications with high repetition rates, pulse picking and regenerative amplifiers
- Two crystals in compensation layout
- Wavefront Distortion: $< \lambda/4$
- Damage threshold: $> 600 \text{ MW/cm}^2$ at 1064 nm, 10 ns, 1 Hz (typical, not guaranteed)
- SC version with short crystals
- Optionally available with integrated Brewster polarizer: RTPC n BP
- Optionally available with integrated $\lambda/4$ disk: RTPC n WP
- All order numbers valid for 1064 nm
- Other specifications on request
- Please state the applied wavelength when ordering.

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
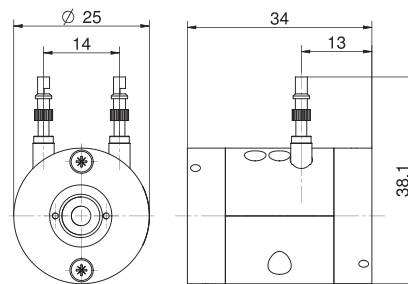
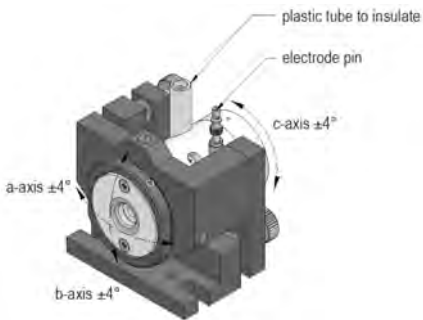
Pockels Cells Positioner


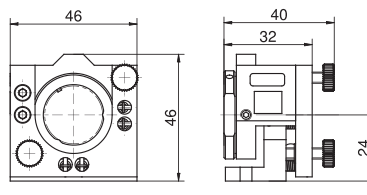
- Compact and stable design
- Easy adjustment of pitch, yaw and rotation
- Adjustment via fine-threaded screws
- For pockels cells with a diameter up to 35 mm
- Optional special OEM modifications available

Positioner 25

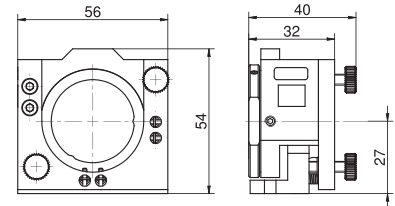Positioner 35

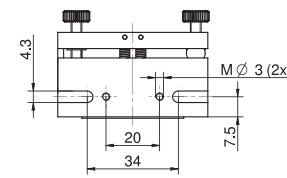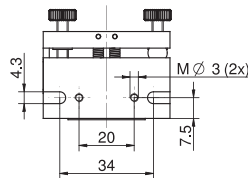Pockelszellen-Positionierer Pockels Cells Positioner

Produktbezeichnung	Abmessungen (mm)	Strahlhöhe (mm)	Bereich Verkippung	Durchmesser Pockelszelle (mm)	Part No.
Item Title	Dimensions (mm)	Beam height (mm)	Tilt range	Diameter pockels cell (mm)	Part No.
Positioner 25	46 x 46 x 40	24	±4	12.7	84503021127
Positioner 25	46 x 46 x 40	24	±4	19	84503021190
Positioner 25	46 x 46 x 40	24	±4	21	84503021210
Positioner 25	46 x 46 x 40	24	±4	23	84503021230
Positioner 25	46 x 46 x 40	24	±4	25	84503021250
Positioner 25	46 x 46 x 40	24	±4	25.4	84503021254
Positioner 35	56 x 54 x 40	24	±4	35	84503021350

Instruments

High-quality laser and light sources for
demanding applications

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Our quality criteria:

- *RoHS and CE-compliant*
- *Laser manufactured in 100 % cleanroom atmosphere*
- *Laser modules sealed in protective atmosphere*
- *Long service life*

LINOS Laser

"A laser is a solution seeking a problem" (Theodore Maiman)

The world's first working laser was invented by U.S. physicist Theodore Maiman in 1960. Today, there are countless applications that simply could not exist without laser technology. Qioptiq helps to make it all possible, with a broad spectrum of laser types for wavelengths ranging from 375 nm to 1064 nm, and output levels of up to 800 mW.

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Special Features:

Our HeNe laser Series can be integrated in our mounting systems directly or with adapters.

Ideal Areas:

Microscopy; endoscopy; quality assurance; inspection systems; image processing; life sciences; biology; Metrology; spectroscopy; medical technology; bioanalysis; fluorescence excitation; materials testing and processing; sensor technology; telecommunications.

Inhalt

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iFLEX-iRIS Laser	562
iFLEX-iRIS Fiber Coupled Laser	564
iFLEX-iRIS CDRH Compliant Interlock	566
iFLEX-iRIS OEM power supply	567
iFLEX-iRIS Accessories	567

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Content

Diode and DPSS Laser

<i>iFLEX-iRIS Laser</i>	562
<i>iFLEX-iRIS Fiber Coupled Laser</i>	564
<i>iFLEX-iRIS CDRH Compliant Interlock</i>	566
<i>iFLEX-iRIS OEM power supply</i>	567
<i>iFLEX-iRIS Accessories</i>	567

HeNe Laser and Accessories

<i>Set Laser and Power Supply: Cylindrical HeNe Lasers with Shutter, Outer Diameter 35 and 45 mm</i>	568
<i>Set laser and power supply: Cylindrical HeNe Lasers, Outer Diameter 31.7 and 44.25 mm</i>	570
<i>Cylindrical HeNe Lasers w/o Power Supply, with Shutter, Outer Diameter 35 and 45 mm</i>	571
<i>Cylindrical HeNe Lasers w/o Power Supply, Outer Diameter 31.7 and 44.25 mm</i>	572
<i>Power Supply for Cylindrical HeNe Lasers</i>	573
<i>C-Mount Adapter for HeNe-Laser with Outer Diameter of 35 mm</i>	574
<i>C-Mount Adapter for HeNe Lasers with Outer Diameter of 45 mm</i>	574
<i>Laser Adapter C 35 for HeNe Laser with Outer Diameter 31.7 mm</i>	575
<i>Microbench Adapter for HeNe-Laser with 31.7 mm Outer Diameter</i>	575
<i>Beam Blocking Shutter, Mechanical</i>	576
<i>Adapter FIC / Microbench</i>	576

iFLEX Agile cw OPO

<i>iFLEX-Agile Series cw-OPO</i>	577
----------------------------------	-----

iFLEX-iRIS with £1 coin

iFLEX-iRIS Laser

The iFLEX-iRIS is a high performance solid-state laser system. All the smart control electronics and thermoelectric cooling (TEC) are fitted inside this small laser. The design includes automatic closed loop control in cw mode, for excellent long term power stability. Ultra-stable performance is also achieved when the laser is in modulation mode. This allows you to achieve measurements that are precise and repeatable, day after day.

Easy to use:

All wavelengths in the same compact size with the same control interface. Performance Guarantee: All lasers are manufactured in clean-rooms, in our ISO9001 compliant factory. Each laser is tested before dispatch, and supplied with a certificate of conformance. Qioptiq design and build lasers under the guidelines given in the international telecommunication standard, Telcordia GR468. This high test standard and ongoing reliability program gives our customers confidence in our product reliability.

Key Features:

- Same compact design for all wavelength options
- Integrated control electronics
- Exceptional power stability
- Low amplitude noise
- Output beam 0.7 ± 0.2 mm
- Operating mode cw

Options:

- CW with output power adjustment options
- Modulation options
- Fiber coupled options

CDRH compliant:

A power supply is required for operation and is sold separately. The iFLEX-iRIS CDRH interlock power supply is required for CDRH compliant systems; and it is also recommended for general laboratory use. An OEM power supply is also available, or the iFLEX-iRIS may be powered and controlled directly through an instrument's power bus.

Applications include:

Metrology, flow cytometry, confocal and laser scanning microscopy, dynamic light scattering, forensics, TIRF, high throughput screening, high-resolution imaging systems, DNA screening, genomics, ophthalmology and more.

Accessories:

- CDRH remote interlock with power supply
- OEM power supply
- kineMATIX-P2 KIT (required for coupling kine FLEX fiber optic to iFLEX-iRIS)
- kineFLEX single-mode polarization maintaining fiber (see kineFLEX)
- Contact us for pricing on other wavelengths, power levels, modulation and fiber options. Standard lead time: 8 weeks.

iFLEX-iRIS Laser

<i>Item Title</i>	<i>Wavelength (nm)</i>	<i>Output power (mW)</i>	<i>Power adjustment (%)</i>
iFLEX-iRIS-X0-375-0.7-50-NP	375	50	Off, 0.1-100%
iFLEX-iRIS-X0-405-0.7-50-NP	405	50	Off, 0.1-100%
iFLEX-iRIS-X0-445-0.7-75-NP	445	75	Off, 0.1-100%
iFLEX-iRIS-X0-458-0.7-70-NP	458	70	Off, 0.1-100%
iFLEX-iRIS-X0-473-0.7-75-NP	473	75	Off, 0.1-100%
iFLEX-iRIS-X0-488-0.7-40-NP	488	40	Off, 0.1-100%
iFLEX-iRIS-X0-505-0.7-50-NP	505	50	Off, 0.1-100%
iFLEX-iRIS-X0-515-0.7-50-NP	515	50	Off, 0.1-100%
iFLEX-iRIS-X0-520-0.7-30-NP	520	30	Off, 0.1-100%
iFLEX-iRIS-X0-532-0.7-40-NP	532	40	Off, 50-100%
iFLEX-iRIS-X0-561-0.7-40-NP	561	40	Off, 15-100%
iFLEX-iRIS-X0-594-0.7-20-NP	594	20	Off, 50-100%
iFLEX-iRIS-X0-633-0.7-70-NP	633	70	Off, 0.1-100%
iFLEX-iRIS-X0-637-0.7-100-NP	637	100	Off, 0.1-100%
iFLEX-iRIS-X0-642-0.7-100-NP	642	100	Off, 0.1-100%
iFLEX-iRIS-X0-660-0.7-80-NP	660	80	Off, 0.1-100%
iFLEX-iRIS-X0-670-0.7-10-NP	670	10	Off, 0.1-100%
iFLEX-iRIS-X0-730-0.7-20-NP	730	20	Off, 0.1-100%
iFLEX-iRIS-X0-780-0.7-70-NP	780	70	Off, 0.1-100%
iFLEX-iRIS-X0-830-0.7-100-NP	830	100	Off, 0.1-100%
iFLEX-iRIS-X0-852-0.7-35-NP	853	35	Off, 0.1-100%

iFLEX-iRIS fiber coupled laser using kineFLEX

iFLEX-iRIS Fiber Coupled Laser

The iFLEX-iRIS is a high performance solid-state laser system. All the smart control electronics and thermoelectric cooling (TEC) are fitted inside this small laser. The design includes automatic closed loop control in cw mode, for excellent long term power stability. Ultra-stable performance is also achieved when the laser is in Modulation mode. This allows you to achieve measurements that are precise and repeatable, day after day.

Fiber coupled output:

A kineFLEX fiber delivery system is included with the fiber coupled versions of the iFLEX-iRIS laser. This single-mode polarization maintaining fiber optic output reduces beam wander to $<1\mu\text{rad}/\text{oC}$ and provides a spatially filtered, near-perfect Gaussian output beam. Collimated and connector output options are available.

Easy to use:

All wavelengths in the same compact size with the same control interface.

Performance guarantee:

All lasers are manufactured in cleanrooms, in our ISO9001 compliant factory. Each laser is tested before dispatch, and supplied with a certificate of conformance. Qioptiq design and build lasers under the guidelines given in the international telecommunication standard, Telcordia GR468. This high test standard and ongoing reliability program gives our customers confidence in our product reliability.

Key Features:

- Same compact design for all wavelength options
- Integrated control electronics
- Exceptional power stability
- Low amplitude noise
- Operating mode cw

Options:

- CW with output power adjustment options
- Modulation options
- Free space (no fiber) options

CDRH compliant:

A power supply is required for operation and is sold separately. The iFLEX-iRIS CDRH interlock power supply is required for CDRH compliant systems; and it is also recommended for general laboratory use. An OEM power supply is also available, or the iFLEX-iRIS may be powered and controlled directly through an instrument's power bus.

Applications include:

metrology, flow cytometry, confocal and laser scanning microscopy, dynamic light scattering, forensics, TIRF, high throughput screening, high-resolution imaging systems, DNA screening, genomics, ophthalmology and more.

Accessories:

- CDRH remote interlock with power supply
- OEM power supply
- kineMATIX-P2 KIT FOR iFLEX-iRIS
- kineFLEX single-mode polarization maintaining fiber (see kineFLEX)

- **Contact us for pricing on other wavelengths, power levels, modulation and fiber options. Standard lead time: 8 weeks.**

iFLEX-iRIS Laser with 2m SM/PM Fiber with 0.7mm diameter collimated output

<i>Item Title</i>	<i>Wavelength (nm)</i>	<i>Output power (mW)</i>
iFLEX-iRIS-P-2-375-0.7-30-NP	375	30
iFLEX-iRIS-P-2-405-0.7-30-NP	405	30
iFLEX-iRIS-P-2-445-0.7-50-NP	445	50
iFLEX-iRIS-P-2-488-0.7-20-NP	488	20
iFLEX-iRIS-P-2-515-0.7-40-NP	515	40
iFLEX-iRIS-P-2-532-0.7-25-NP	532	25
iFLEX-iRIS-P-2-561-0.7-30-NP	561	30
iFLEX-iRIS-P-2-642-0.7-65-NP	642	65

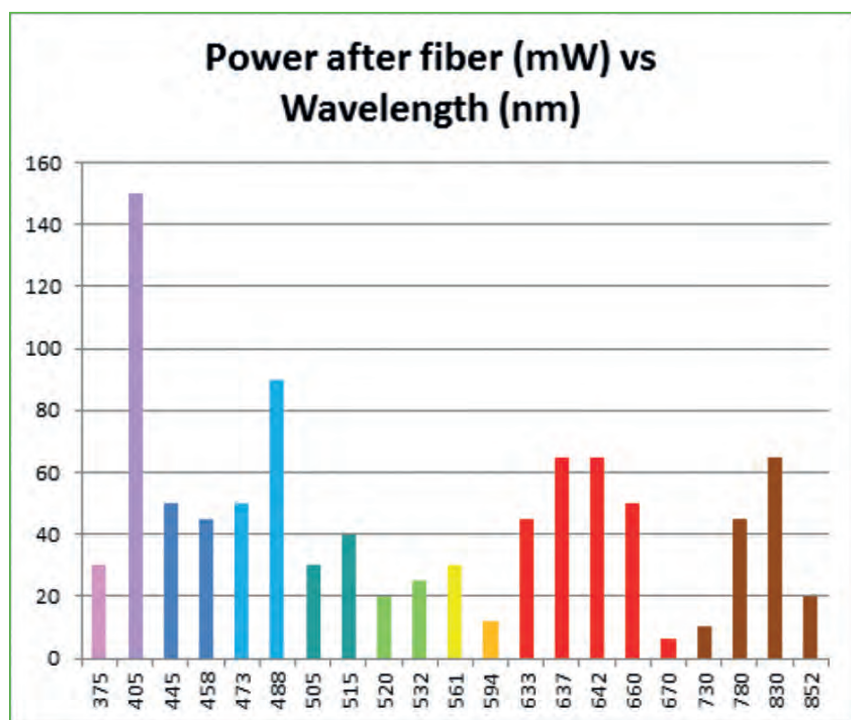
Photo shows 2 pcs. iFLEX-iRIS CDRH compliant interlocks and power supplies

iFLEX-iRIS CDRH Compliant Interlock


The iFLEX-iRIS CDRH compliant interlock remote with power supply unit is recommended for general laboratory use, and complies with CDRH regulations.

It includes a laser safety keyswitch (with 2 pcs. keys) and a laser safety interlock plug, plus LED indicator to show laser emission status. It is provided with a power supply.

Please note that a standard IEC power cable must be acquired separately. This can easily be purchased locally to match your local power plug requirements.

iFLEX-iRIS CDRH Compliant interlock

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
iFLEX-iRIS CDRH compl. interlock	on request

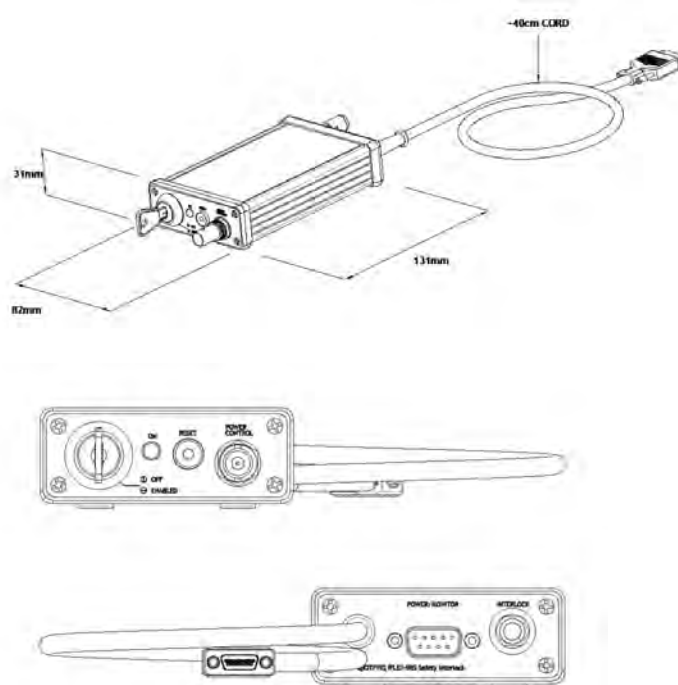iFLEX-iRIS OEM power supply

Qioptiq is pleased to offer this small OEM power supply for the iFLEX-iRIS laser systems.


iFLEX-iRIS lasers deliver optimum performance when they are connected to a clean power supply. For this reason Qioptiq offer either the iFLEX-iRIS OEM or iFLEX-iRIS CDRH compliant power supplies.

Please be aware that this OEM power supply does not include laser safety features, and it is not compliant to CDRH regulations for general laboratory use. Please note that a standard IEC power cable must be acquired separately. This can easily be purchased locally to match your local power plug requirements.

When integrated into OEM systems then power may be supplied to iFLEX-iRIS lasers directly by the instrument.

iFLEX-iRIS OEM Power Supply

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
iFLEX-iRIS OEM power supply	on request

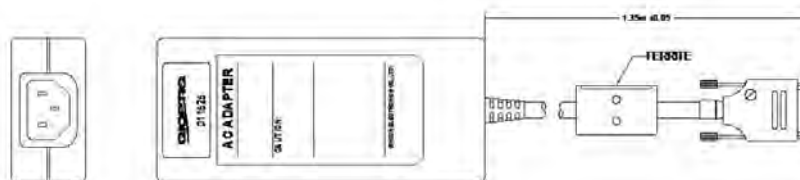iFLEX-iRIS Accessories

Produktbezeichnung	Part No.
<i>Item Title</i>	<i>Part No.</i>
kineMATIX-P2 kit for iFLEX-iRIS	on request
iFLEX-iRIS, optical table mount, 40 mm	on request
iFLEX-iRIS, rail system FLS 40 adapter	on request
iFLEX-iRIS, Microbench adapter	on request
iFLEX-iRIS, Nanobench adapter	on request

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Set Laser and Power Supply: Cylindrical HeNe Lasers with Shutter, Outer Diameter 35 and 45 mm


- Emit visible beam at $\lambda = 543, 594, 633 \text{ nm}$
- Output power from 0.5 up to 20 mW
- Mechanical beam-blocking shutter
- Different versions with polarized laser beam (P)
- Singlemode and multimode versions
- High beam-pointing stability even at elevated ambient temperature
- Power supply, switchable $115 \text{ V} / 230 \text{ V}$
- Key switch
- Emission control
- Connectors for external safety loop
- Output power versions from $25-30 \text{ mW}$ available on request.

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Produktbezeichnung	Wellenlänge (nm)	Ausgangsleistung (mW)	Strahl-Ø ($1/e^2$) (mm)	Strahl-divergenz (mrad)	Transversale Modenstruktur (*)	Polarisation
Item Title	Wavelength (nm)	Output power (mW)	Beam-Ø ($1/e^2$) (mm)	Beam divergence (mrad)	Transverse mode structure (*)	Polarization
Laser Head + Powersupply HeNe 633-0.5 P	632.8	0.5	0.49	1.75	SM	>500:1
Laser Head + Powersupply HeNe 633-0.6	632.8	0.6	0.49	1.7	SM	N/A
Laser Head + Powersupply HeNe 633-1	632.8	1	0.49	1.7	SM	N/A
Laser Head + Powersupply HeNe 633-2	632.8	2	0.75	1.2	SM	N/A
Laser Head + Powersupply HeNe 633-2 P	632.8	2	0.75	1.2	SM	>500:1
Laser Head + Powersupply HeNe 633-5	632.8	5	0.8	1.1	SM	N/A
Laser Head + Powersupply HeNe 633-5 P	632.8	5	0.8	1.1	SM	>500:1
Laser Head + Powersupply HeNe 633-5 M	632.8	5	1.4	4.5	MM	N/A
Laser Head + Powersupply HeNe 633-10	632.8	10	0.7	1.4	SM	N/A
Laser Head + Powersupply HeNe 633-10 P	632.8	10	0.7	1.4	SM	>500:1
Laser Head + Powersupply HeNe 633-10 M	632.8	10	1.5	4	MM	N/A
Laser Head + Powersupply HeNe 633-15	632.8	15	1	1	SM	N/A
Laser Head + Powersupply HeNe 633-15 P	632.8	15	1	1	SM	>500:1
Laser Head + Powersupply HeNe 633-18	632.8	18	1	1	SM	N/A
Laser Head + Powersupply HeNe 633-20	632.8	20	1	1	SM	N/A
Laser Head + Powersupply HeNe 633-18 P	632.8	18	1	1	SM	>500:1
Laser Head + Powersupply HeNe 543-0.5 P	543.5	0.5	0.9	1	SM	>500:1
Laser Head + Powersupply HeNe 543-0.75 P	543.5	0.75	0.85	0.9	SM	>500:1
Laser Head + Powersupply HeNe 543-1	543.5	1	0.88	0.85	SM	N/A
Laser Head + Powersupply HeNe 543-1 P	543.5	1	0.85	0.91	SM	>500:1
Laser Head + Powersupply HeNe 543-1.2 P	543.5	1.2	0.85	0.91	SM	>500:1
Laser Head + Powersupply HeNe 543-1.5	543.5	1.5	0.88	0.85	SM	N/A
Laser Head + Powersupply HeNe 543-2	543.5	2	0.88	0.85	SM	N/A
Laser Head + Powersupply HeNe 543-2.5	543.5	2.5	0.88	0.85	SM	N/A
Laser Head + Powersupply HeNe 594-2 P	594.1	2	0.79	0.95	SM	>500:1

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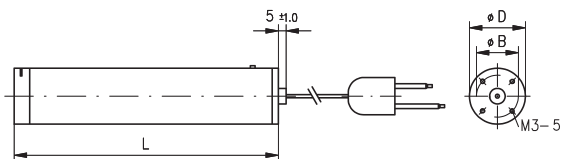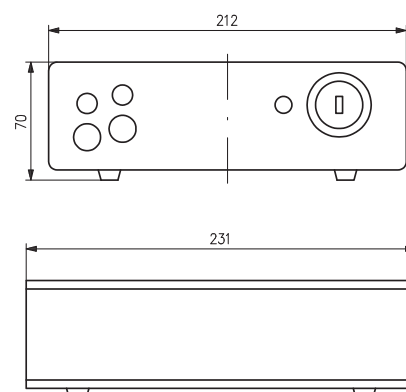
- Transversale Modenstruktur: Singlemode

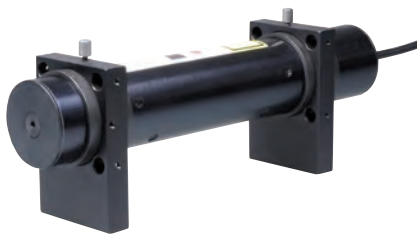
- Schutzklasse IEC Europe: 3R
- Schutzklasse CDRH USA: IIIa

Set laser and power supply: Cylindrical HeNe Lasers, Outer Diameter 31.7 and 44.25 mm

- Emit visible beam at $\lambda=543$ or 633 nm
- Output power from 0.5 up to 2 mW
- Version 543 nm with mechanical beam-blocking shutter
- Different versions with polarized laser beam (P)
- High beam-pointing stability even at elevated ambient temperature
- Power supply, switchable 115 V / 230 V
- HeNe-Laser with outer diameter 31.7 mm fits to LINOS Microbench with adapter G040 691 000
- Power supply, switchable 115 V / 230 V
- Key switch
- Emission control
- Connectors for external safety loop


- Transverse mode structure: Single mode

- Safety class IEC Europe: 3R
- Safety class CDRH USA: IIIa

HeNe-Laser with outer diameter of 31.7 mm attached in LINOS Microbench

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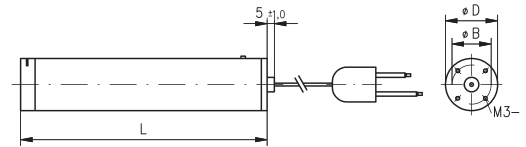Cylindrical HeNe Lasers w/o Power Supply, with Shutter, Outer Diameter 35 and 45 mm

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- Emit visible beam at $\lambda = 543, 594, 633$ nm
- Output power from 0.5 up to 20 mW
- Mechanical beam-blocking shutter
- Different versions with polarized laser beam (P)
- Singlemode and multimode versions
- High beam-pointing stability even at elevated ambient temperature
- Separate power supply, switchable 115 V / 230 V (not included)
- Output power versions from 25-30 mW available on request.

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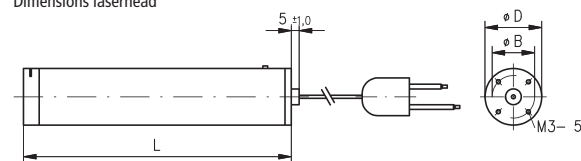
Cylindrical HeNe Lasers w/o Power Supply, Outer Diameter 31.7 and 44.25 mm

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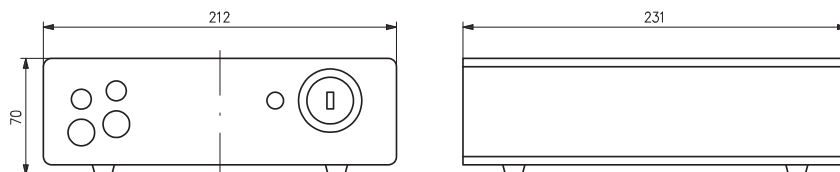
Power Supply for Cylindrical HeNe Lasers


- Key switch
- Emission control
- Connectors for external safety loop
- Input voltage (optional) 230 / 115 VAC

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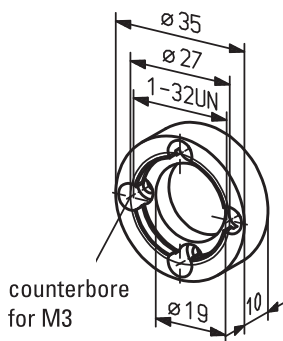
C-Mount Adapter for HeNe-Laser with Outer Diameter of 35 mm

- For assembly of components with a C-Mount thread like beam expanders to a cylindrical HeNe-Laser with 35 mm outer diameter

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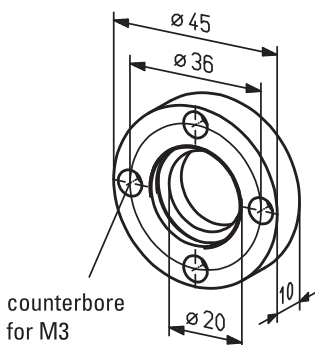C-Mount Adapter for HeNe Lasers with Outer Diameter of 45 mm

- For assembly of components with a C-Mount thread like beam expanders to a Cylindrical HeNe-laser with 45 mm outer diameter

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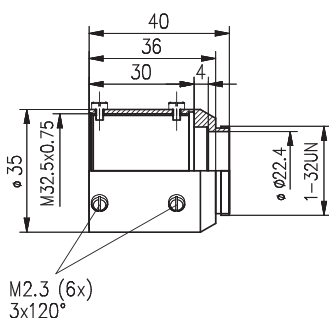
Laser adapter C 35

Laser adapter C 35 with screw-on C-Mount Microbench adapter

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
Laser Adapter C 35 for HeNe Laser with Outer Diameter 31.7 mm

- To mount components with C-Mount thread (e.g. Laser expander systems) to cylindrical HeNe Laser with 31.7 mm outer diameter

Laseradapter C 35 Laser Adapter C 35


Produktbezeichnung	Part No.
Item Title	Part No.
Laser adapter C 35	G040102000

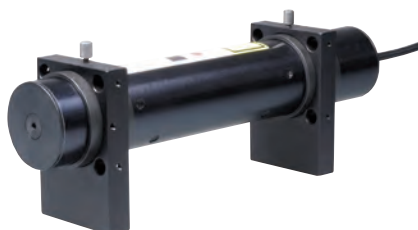HeNe-Laser with outer diameter of 31.7 mm attached in Microbench

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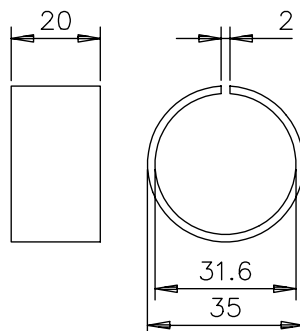Microbench Adapter for HeNe-Laser with 31.7 mm Outer Diameter

- Adapter to attach a cylindrical HeNe Laser with 31.7 mm outer diameter to Microbench Mounting plate 35 (G06 1047 000) and Holder 35 (G06 5084 000).

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
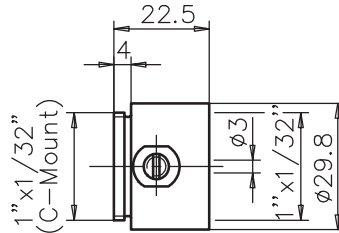
Beam Blocking Shutter, Mechanical

- C-Mount thread
- Direct connection to tube system C possible
- Reduction adapter available for fitting to Microbench

Strahlunterbrecher, mechanisch (Shutter) Beam blocking shutter, mechanical

Produktbezeichnung	Part No.
Item Title	Part No.
Beam blocking shutter, mechanical	G040212000

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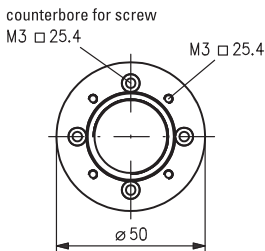Adapter FIC / Microbench

- 30 mm flange diameter for installing in e.g. LINOS Microbench mounting plate 30 or holder 30
- With 4x M3 threads directly attachable to mounting bracket DPL, G04 0872 000
- With internal C-Mount thread

Adapter FIC / Mikrobank Adapter FIC / Microbench


Produktbezeichnung	Part No.
Item Title	Part No.
Adapter FIC/Microbench	G173013000

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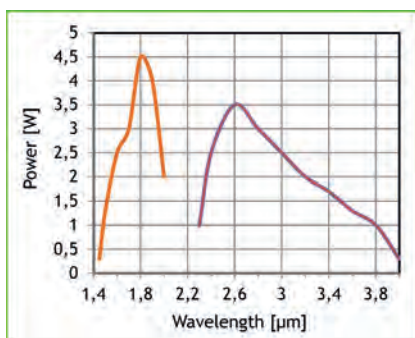iFLEX-Agile Series cw-OPO

The New iFLEX-Agile Series comprises various combinations of continuous-wave optical parametric oscillator (cw-OPO) modules and pump-sources featuring exceptionally broad wavelength tunability.

The modular concept allows modification of wavelength tuning ranges, emission linewidths and output power levels to cost-efficiently meet the custom requirements for specific applications.

Key Features

- Broad wavelength tunability at high-power continuous-wave outputs >1W
- Rapid tunability via USB interface across 1480-2000nm and 2300-3800nm from a single OPO-module without exchange of optics
- Exceptional emission beam quality

Options

- Narrow linewidths <1MHz over the whole emission range
- Higher power levels
- Broader wavelength emission range, e.g. additional 740-950nm
- Wavelength tuning options

Applications

- Trace-gas analysis
- Spectroscopy
- Metrology
- Detector calibration
- Mass spectrometry
- Material (specification) tests

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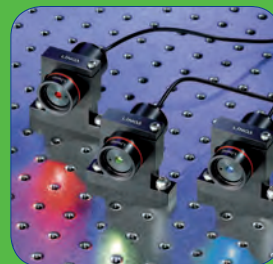
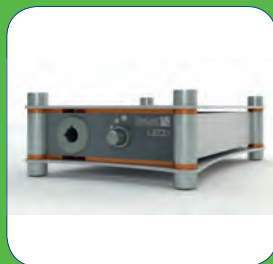
- Made in Germany
- RoHS- und CE-konform
- Lange Lebensdauer

Our quality criteria:

- *High-quality German engineering*
- *RoHS and CE-compliant*
- *Long service life*

LINOS Light Sources

Qioptiq provides a wide variety of light sources. We offer the classic broadband light sources in addition to monochromatic laser light sources. Moreover, Qioptiq has halogen and xenon fiber light sources and LED light sources for reproduction with a CCD camera or observation unit. Our small halogen lamp housings, directional lamps and spectral lamps – compatible with all our assembly components – are especially well suited for laboratory setups.

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Special Features:

We have a broad range of accessories such as light guides, focusing methods and filters for all cold light sources. Condensers, fans and power supplies complete the line of lamp housing products.

Ideal Areas:

Microscopy; endoscopy; quality assurance; inspection systems; image processing; life sciences; biology

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Content

Microbench Light Sources	
<i>Tungsten-Halogen Lamp House, 50 W</i>	582
<i>Tungsten-Halogen Lamp Socket</i>	582
<i>Tungsten-Halogen Lamp 12 V</i>	582
<i>LED Light Sources HighLED</i>	583
<i>Condenser C35 with Ancillary Lens</i>	584
<i>Dispersion Plate in Retaining Mount</i>	584
<i>Polarization Filter in Retaining Mount</i>	585
Lamphouse IL and Accessories	
<i>Lamphouse IL</i>	586
<i>Spectral Lamps</i>	587
<i>Mercury Arc Lamp</i>	588
<i>Xenon Arc Lamp</i>	589
<i>Halogen Lamp Power Supply PSH 100</i>	590
<i>Spectral Lamp Power Supply PSP 60</i>	591
<i>Mercury Arc Lamp Power Supply PSQ 50AC</i>	592
<i>Xenon Lamp Power Supply PSX 75</i>	593
<i>Mounting Bracket 40/65</i>	594
<i>Microbench Adapter IL</i>	594
<i>Condenser; Two-Lens; 105-385 mm; VIS</i>	595
<i>Condenser; Two-Lens; 105-385 mm; UV</i>	595
<i>Condenser; One-Lens 220 mm to ∞; VIS</i>	596
<i>Condenser; One-Lens 220 mm to ∞; UV</i>	596
Fiber Illumination System	
<i>Fiber Illumination Systems, Overview</i>	597
<i>Flexible Fiber Light Guide</i>	598
<i>Focusing Lens</i>	598
<i>Fiber Optic Ring Light</i>	599
<i>Focusing lens</i>	600
<i>Liquid Fiber Light Guides LLG VIS</i>	600
<i>Liquid Fiber Light Guides LLG IR</i>	601
<i>Liquid Fiber Light Guides LLG UV</i>	602
<i>Focusing Lens Attachments FA-LLG</i>	603
LED Light Sources	
<i>Cold Light Sources CLS-LED3</i>	604
<i>Cold Light Source CLS-LED5</i>	605
<i>Cold Light Sources CLS-LED7</i>	606
<i>LED-Industry Lamp IL100-c</i>	607
<i>Modules for IL100-c</i>	608
<i>Microbench Light Source LED ML</i>	609
X-Cite® Fluorescence Illumination Solutions	
<i>X-Cite® Product Overview</i>	610
<i>Fluorescence Microscope Light Sources</i>	610
<i>Power Measurement in Fluorescence Imaging</i>	610
Accessories	
<i>LED coupler</i>	612

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Tungsten-Halogen Lamp House, 50 W

- Uses 12 V / 50 W Tungsten-Halogen Lamp (not included)
- Includes Concave Mirror (G06 03611 000) and Aspheric Condenser (G06 3097 000)
- Ø 4 mm female plugs for connecting electrical supply

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Tungsten-Halogen Lamp Socket


- Component of G06 5005 000 tungsten-halogen lamp house
- Concave mirror and tungsten-halogen lamp not included
- Ø 4 mm female plugs for connecting electrical supply
- Mounting options:
- M4/M6 tapped holes for mounting on threaded posts
- Mounts on breadboards, mounting bases, baseplates
- Mounts on Microbench rods
- Tungsten-halogen lamp 12 V/50W (G065031000), see below.

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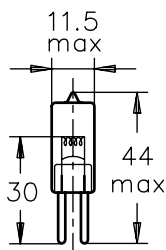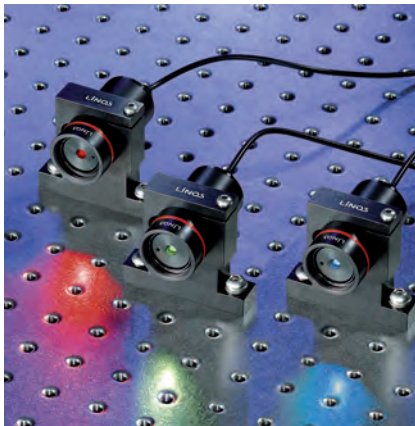
G12 7435 000

HighLED Light Sources in Clamp Holder 30 (G06 1243 000, not included)

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LED Light Sources HighLED



- Super bright LED light sources
- Brightness adjustable
- Current stabilized
- Connection of Qioptiq standard optics with C-Mount
- Compatible to the LINOS Microbench system
- LED: Luxeon Star
- Supply Voltage: 12 V / DC (7 V min, 13 V max)
- Supply Voltage Connection: Binder serie 719
- Current Stabilization: 150 mA ± 2%
- Current Control: 0-150 mA
- Operating Temperature: 0-40 °C
- Mount, Diameter: 30 mm (2.56 in.)
- Mount, Total Length: 65 mm (2.56 in.)
- Weight: 60 g (2.1 oz.)
- A wide range of accessories to the LED light sources HighLED can be found in the following.


LED Light Sources HighLED LED Light Sources HighLED

Produktbezeichnung	Farbe	Typ. Wellenlänge	Typ. Lichtleistung	Artikel-Nr.
Item Title	Color	Typ. wavelength	Typ. output power	Part No.
HighLED-W	white	5550 K	4.0 mW at 440 nm	G065150000
HighLED-B	blue	470 nm	2.5 mW at 470 nm	G065151000
HighLED-G	green	530 nm	2.5 mW at 530 nm	G065152000
HighLED-R	red	630 nm	4.0 mW at 630 nm	G065153000

LED Light Sources HighLED Spannungsversorgung LED Light Sources HighLED power supply

Produktbezeichnung	Input	Output	Artikel-Nr.
Item Title	Input	Output	Part No.
Power supply	100-240 V AC / 50-60 Hz	12 V DC / 450 mA	G127435000

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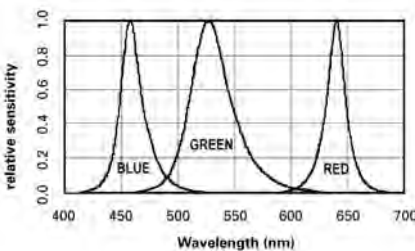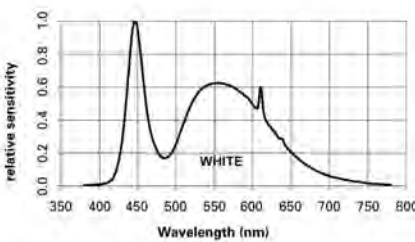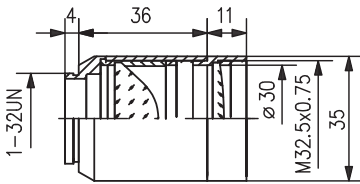
Condenser C35 with Ancillary Lens

HighLED Accessory

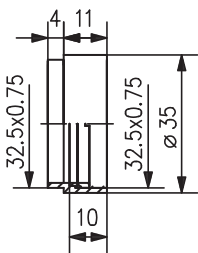- Collimating configuration for HighLED
- Compatible to the LINOS Systems Tube Mounting System C and Microbench
- Easy to detach to assemble other optical configurations
- Thread M32.5x0.75 to attach other ancillary components e.g. polarizers etc.
- Aspheric condenser with plano-convex lens
 $f = 200 \text{ mm}$
- C-Mount thread
- Outer Diameter 35 mm

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
Dispersion Plate in Retaining Mount

HighLED Accessory

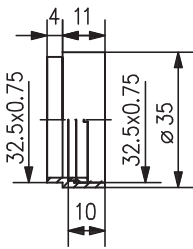- Compatible to the LINOS Systems Tube Mounting System C and Microbench
- Ground glass for diffuse light
- Clear Aperture $\varnothing 30 \text{ mm}$
- Outer Diameter 35 mm

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Polarization Filter in Retaining Mount


HighLED Accessory

- To create linear polarized light
- Compatible to the LINOS Systems Tube Mounting System C and Microbench
- Can be combined with other retaining mounts
- Combine with a second polarizer to vary the light intensity
- Outer Diameter 35 mm

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Polarization filter in retaining mount

Produktbezeichnung	Artikel-Nr.	Preis (Euro) (gültig für DE, AT, CH)
Item Title	Part No.	Price (Euro) (valid in DE, AT, CH)
Polarization filter in retaining mount	G065245901	75,00

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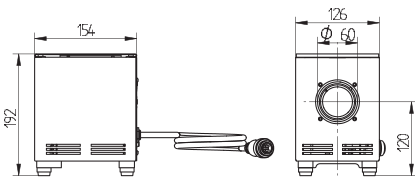Lamphouse IL

The IL Series Lamphouse from Qioptiq consists of a universal lamp housing for halogen, spectral, xenon and mercury vapor lamps. The combination with an electronic lamp power supply from the PS Series gives you a high-quality, sophisticated lighting system for your applications.

The universal lamp housing is exceptionally compact, and the precision adjustment of its mechanical components makes your lighting configurations easy to reproduce. The lamp socket, mirrors and collector can be adjusted from outside the housing, using the adjustments on the back panel. Even when operating under thermal load, the mechanical positioning remains extremely stable. Openings in the top and bottom housing panels allow air to circulate around the lamp without permitting any light to escape. This noiseless and vibration-free convection cooling helps ensure the conditions for stable light emission from the lamp. A $\varnothing 60$ mm collector guide holds the condensers. The lamp housings come with four screw-in rubber feet as standard equipment, any of which can be replaced by an adjustment foot if desired.

- Lamp housings for halogen and spectral lamps, Hg and Xe
- Igniter built in to the lamp housing for Xenon lamps
- External, thermally stable and reproducible mechanical adjustment of lamps, mirrors and collectors
- Easy lamp replacement
- Light-tight
- Splinter shield included in delivery
- Safety-circuit function
- Comprehensive range of accessories
- Compatible with LINOS Microbench

- Housing: Formed sheet metal housing, light-tight, textured paint: white (RAL 9003)
- Dimensions (HxWxD): 169 x 126 x 154 mm (6.65 x 4.96 x 6.06 in.)
- Weight: 3.4 kg (7.5 lb)
- Clearance, Optical Axis – Lower Edge of Lamp Housing : 120 mm (4.72 in.)
- Fastening, Collector: Collector guide with z-adjustment
- Fastening, Lamp: Lamp adapter with X-Y adjustment
- Fastening, Microbench Adapter: M4 hole circle, 73 mm
- Reflector: Mirror with X-Y-Z adjustment
- Lamp Socket: Lamp adapter (lamp-type specific)
- Power Supply: Electronic ballast (lamp specific)
- Electrical Connection: Cable, 1.5 m long; electrical connection permanently attached to lamp housing, plug coded to match electronic ballast
- Ambient Temperature: 0 °C to + 40 °C
- Storage Temperature: - 40 °C to +70 °C
- Device Protection Class: I
- Type of Protection: IP 20 (EN 60529)
- Safety Circuit: Shut off when device is opened

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Spectral Lamps

Spectral lamps are discharge lamps that emit the line spectrum of inert gases and metal vapors with high luminous density or radiance. They are utilized wherever a line spectrum or monochromatic radiation is required; primarily in the fields of optics, radiation physics, spectroscopy, chemistry and medicine.

- *Nominal current: 1 A*
- *Socket: Pico 9*
- *Burning position: Vertical*

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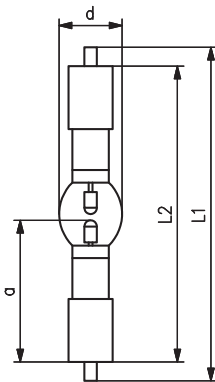
Mercury Arc Lamp


HBO lamps are short-arc lamps with a discharge arc that burns in a mercury vapor atmosphere under high pressure. Typical areas of application include fluorescence microscopy, fluorescence endoscopy, schlieren photography and hologram reduction.

The HBO 50 W/AC is solely an alternating-current lamp and must not be connected to direct current.

- Ideal lamp for Lamphouse ILQ 50AC for Mercury Arc Lamps G03 0532 000
- High radiance
- High optical quality
- Multiple-line spectrum
- High radiation output in the UV and visible range
- Long service life
- Power rating: 50 W
- Voltage drop: 39-45 V
- Nominal current: 1.3 A
- Luminous flux: min. 1650 lm
- Luminosity: min. 200 cd
- Luminous density: min. 27000 cd/cm²
- Luminous field: 0.3 x 1.0 mm
- Service life: 100 h
- Burning position: vertical to 45 °
- Socket: SFa 6-2

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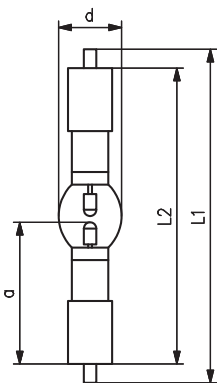Xenon Arc Lamp

XBO xenon lamps are short arc lamps in which the luminous arc burns in an atmosphere of pure xenon gas at high pressure. XBO lamps are used in scientific and technical applications such as:

- Fluorescence Spectrometers
- Light sources for microscopy / fluorescence microscopy
- Light sources for endoscopy / fiber optics
- Graphic scanners
- Color matching lights
- Color measuring units

- Ideal lamp for Lamphouse ILX 75 for Xenon Lamps G03 0534 000
- High luminance
- Daylight color temperature of approx. 6000 K
- Continuous spectrum in the visible range
- Constant color properties throughout the life time of the lamp
- High arc stability
- Almost full luminous flux immediately after ignition
- High color rendering index (Ra > 95)

- Power rating: 75 W
- Voltage drop: 14 V
- Nominal current: 5.4 A
- Luminous flux: 1000 lm
- Luminosity: min. 100 cd
- Luminous density: 40000 cd/cm²
- Luminous field: 0.3 x 0.8 mm
- Service life: 400 h
- Burning position: vertical to 100°
- Socket: Anode SFa 9-2, Cathode SFa 7.5-2

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Halogen Lamp Power Supply PSH 100

- Equipment supplied: power supply with power cord and operating instructions

The PSH 100 Power Supply is intended solely for operation of halogen lamps with max. 12 V/100 W in suitable lamp housings.

The output voltage from the ballast can be adjusted by the user with high precision using the 10-turn potentiometer on the front panel of the device. The potentiometer adjustment knob can be secured to prevent inadvertent changes in the output voltage. The display on the front panel of the instrument can show output voltage or output current, depending on user selection. A service-hour counter is provided to record the number of hours of operation. The counter can be reset to 0 by the user.

- Dimensions (WxHxD): 175 x 115 x 290 mm (6.89 x 4.53 x 11.42 in.)
- Weight: 2.8 kg (6.2 lb)
- Output Voltage; adjustable: 1 to 12.1 VDC
- Line Regulation: $< \pm 0.2 \%$
- Output Voltage Ripple: $< 1 \%$ peak to peak
- Output Current: Max. 10 A
- Short-Circuit Strength: Continuous short-circuit protected
- Input Voltage: 100 to 240 VAC ($\pm 10\%$)
- Frequency: 50 to 60 Hz
- Power Consumption: Max. 150 VA
- Fuse: 2 T4 A; breaking capacity H
- Permissible Ambient Temperature Range (storage): -20 to $85 \text{ }^\circ\text{C}$
- Permissible Relative Humidity (storage): 10 to 90%, non-condensing
- Permissible Ambient Temperature Range (operation): -0 to $40 \text{ }^\circ\text{C}$
- Permissible Relative Humidity (operation): 10 to 90%, non-condensing
- Device Protection Class: I
- Type of Protection: IP 20 (EN 60529)
- Overheating Protection: Overheating causes the equipment to switch off. Once the equipment has cooled down, lamp output switches back on automatically.
- Overload Protection: Output is protected from continuous short-circuit current
- Limitation of Emissions: EN 61326 (Class B)
- Defined Immunity to Interference: EN 61326 Appendix A
- Voltage Fluctuation, Flicker: EN 61000-3-3
- Harmonic Current Emissions: EN 61000-3-2

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Spetral Lamp Power Supply PSP 60

The Power Supply PSP 60 is solely designed to operate the Qioptiq Lamp house ILP 60 (G03 0531 000) using the types of spectral lamps from OSRAM listed below

Approved types of lamps:

- Hg100
- HgCd/10
- He/10
- Na/10
- Ne/10
- Tl/10
- Zn/10
- Cs/10*
- K/10*
- Rb/10*

* Not available by Qioptiq.

- Dimensions (WxHxD): 165 x 85 x 240 mm (6.5 x 3.35 x 9.45 in.)
- Weight: 1,4 kg (3.1 lb)
- Output Current (Nominal value): 1.0 A ± 0.05 A
- Stability (8 h): < 10⁻² Max.
- Output Current (Burn-In Current): max. 1.8 A
- Output Current Pulsation: < 5%
- Burning Voltage (Nominal value): 10 ... 60 VAC
- Open Circuit Voltage in the Ignition Phase (approx. 5 sec.): max. 600 VAC = Ignition Voltage
- Open Circuit Voltage in the Operating Phase: max. 70 VAC (Shutdown)
- Input Voltage: 220 ... 240 VAC (±10%)
- Frequency: 50 ... 60 Hz
- Power Consumption: max. 100 VA
- Fuse: 2x T1,6 A; Breaking Capacity H
- Permissible Ambient Temperature Range (operation): 0 to 40°C
- Permissible Relative Humidity (operation): 10 to 90 %, non-condensing
- Device Protection Class: I
- Type of Protection: IP20 (EN 60529)

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Mercury Arc Lamp Power Supply PSQ 50AC

This electronic AC Power Supply replaces the older-model electrical ballasts and offers outstanding properties which in turn extend service life of the lamp. Moreover, it provides significantly improved light quality, and can be used anywhere in the world with no switching needed. In addition to the traditional areas of use, this considerably smaller and lighter instrument opens up many new possibilities as well, such as in connection with camera applications, for example. The electrical ballast is based on a switching power supply. In comparison to the power supplies with choke coils usually implemented for mercury vapor lamps, this model enables the following advantages:

- *Energy-saving operation enables a lamp service life of approx. 500 h*
- *Greater efficiency*
- *Extremely low ripple of the luminous flux*
- *Wide range input*
- *Economies of weight and volume*

The power supply also generates the starting voltage required to ignite the connected lamp. No additional igniter is needed. The starting voltage is generated automatically when the power supply is switched on, after the safety circuit is closed, and switches off again automatically once the connected mercury short-arc lamp as been successfully ignited.

- *Dimensions (WxHxD): 165 x 86 x 235 mm (6.50 x 3.39 x 9.25 in.)*
- *Weight: 1.4 kg (3.09 lb)*
- *Output Power: 50 ± 3 W*
- *Stability (8 h): ≤ 2 %*
- *Combustion Current: Max. 2.0 A*
- *Combustion Voltage: 33 to 55 V AC*
- *Output Frequency: 151 Hz*
- *Pulse Wave Shape: Rectangular*
- *Luminous Flux Modulation: < 5 %*
- *Ignition: Automatic, with ignition time limit*
- *Ignition Voltage: Approx. 500 VAC eff.*
- *Input Voltage: 100 to 240 VAC (±10%)*
- *Frequency: 50 to 60 Hz*
- *Power Consumption: Max. 150 VA*
- *Fuse: 2 T4 A; breaking capacity H*
- *Permissible Ambient Temperature Range (storage): -20 to 85 °C*
- *Permissible Relative Humidity (storage): 10 to 90%, non-condensing*
- *Permissible Ambient Temperature Range (operation): 0 to 40 °C*
- *Permissible Relative Humidity (operation): 10 to 90%, non-condensing*
- *Device Protection Class: I*
- *Type of Protection: IP 20 (EN 60529)*
- *Overheating Protection: Combustion voltage is cut; output remains connected to the mains*
- *Safety Circuit: Present*
- *Outputs: Connected to the mains*
- *Limitation of Emissions: EN 61326 (Class B)*
- *Defined Immunity to Interference: EN 61326 Appendix A*
- *Voltage Fluctuation, Flicker: EN 61000-3-3*
- *Harmonic Current Emissions: EN 61000-3-2*

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
Xenon Lamp Power Supply PSX 75

The PSX 75 is Power Supply made solely for operation of type XBO 75 W/2 model Xenon short-arc lamps from OSRAM and equivalent types in suitable lamp housings. The corresponding Qioptiq lamp housing, ILX 75, has a built-in igniter that generates the high voltage necessary to ignite the Xenon short-arc lamp. The PSX 75 power supply provides a no-load voltage (UL = 55 to 65 V DC) to the lamp output for operation of the igniter.

- Dimensions (WxHxD): 167 x 85 x 222 mm (6.57 x 3.35 x 8.74 in.)
- Weight: 2.2 kg (4.85 lb)
- Output Power: 86 W
- Output Current: Nominal value 5.4 A ± 0.3 A
- Stability (8 h): ≤ 10-2
- Output Current Ripple (50 Hz to 200 kHz): < 50 mApp
- No-Load Voltage: 55 to 65 VDC
- Combustion Voltage: 10 to 18 VDC
- Short-Circuit Strength: Continuous short-circuit protected
- Input Voltage: 100 to 240 VAC (±10%)
- Frequency: 50 to 60 Hz
- Power Consumption: Max. 130 VA
- Fuse: 2 T2 A; 5 x 20 mm; breaking capacity H
- Permissible Ambient Temperature Range (storage): -20 to 85 °C
- Permissible Relative Humidity (storage): 10 to 90%, non-condensing
- Permissible Ambient Temperature Range (operation): 0 to 40 °C
- Permissible Relative Humidity (operation): 10 to 90%, non-condensing
- Device Protection Class: I
- Type of Protection: IP 20 (EN 60529)
- Overheating Protection: Present
- Safety Circuit: Present
- Overload Protection: The lamp output is continuous short-circuit protected
- Output Power Limit: As the lamp's combustion voltage increases over the service life of the lamp, the power limit rises as well. This is why the device limits the maximum output power to 86 W.

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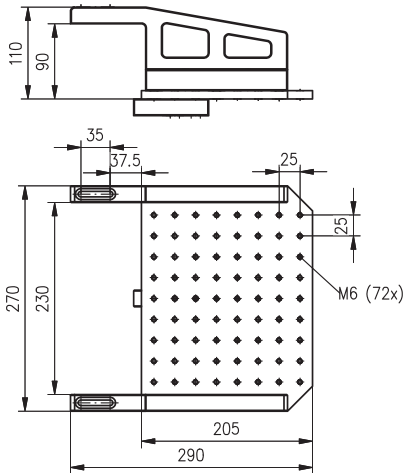
Mounting Bracket 40/65

With the mounting bracket 40/65 the IL series lamp housings can be affixed laterally to an optical table to save valuable space on the work surface. Another practical application is the adaptation of a 120 mm beam height to the standard 40 and 65 mm beam heights of LINOS structural components. This is achieved by simply removing an intermediate piece from the bracket. The base of the mounting bracket has M6 threaded holes (pattern: 25 mm).

- Lateral installation on optical tables
- For adapting beam heights of 120 mm to 40 or 65 mm
- Work surface with M6 threaded holes in a 25 mm pattern

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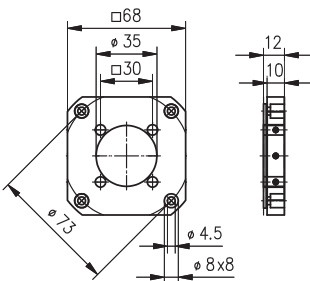Microbench Adapter IL

The Microbench IL adapter is the interface to the Microbench and to Tube Mounting System C from LINOS, placing a broad range of mechanical and optical components at your disposal.

- For easy installation in the LINOS Microbench system
- Free Ø: 35 mm

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Condenser; Two-Lens; 105-385 mm; VIS

- Wavelength range: 420 to 680 nm
- Focal length: 32 mm
- Back focal length: 105 to 385 mm (measured from front panel of lamp housing)
- Numerical aperture: 0.52
- Type: aspherical lens (B270) and plano-convex lens (N-BK7)
- Outer diameter: 60 mm
- Filter thread: M55x0.75

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Condenser; Two-Lens; 105-385 mm; UV

- Wavelength range: 230 to 390 nm
- Focal length: 31 mm
- Back focal length: 105 to 385 mm (measured from front panel of lamp housing)
- Numerical aperture: 0.52
- Type: aspherical lens and plano-convex lens (quartz)
- Outer diameter: 60 mm
- Filter thread: M55x0.75

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Condenser; One-Lens 220 mm to ∞ ; VIS

- Wavelength range: 420 to 680 nm
- Focal length: 40 mm
- Back focal length: 220 mm to infinity
- Numerical aperture: 0.52
- Type: aspherical lens (B270)
- Outer diameter: 60 mm
- Filter thread: M55x0.75

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Condenser; One-Lens 220 mm to ∞ ; UV

- Wavelength range: 230 to 390 nm
- Focal length: 40 mm
- Back focal length: 220 mm to infinity
- Numerical aperture: 0.52
- Type: aspherical lens (fused silica)
- Outer diameter: 60 mm
- Filter thread: M55x0.75

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Fiber Illumination Systems, Overview

The fiber optic cold illuminator consists of a cold light source and a fiber light guide that transmits light from the light source to the object. Our large selection of light guide fiber bundles, liquid light guides with a variety of accessories and the appropriate light sources provide the optimum solutions to lighting problems for the following applications:

- Microscopy
- Endoscopy
- Work place illumination (assembly, precision engineering)
- Quality assurance
- Inspection systems
- Photography
- Adhesive and plastic curing
- Fluorescence excitation
- Irradiation

The essential criteria for selecting a cold light source are the following:

- Intensity output
- Spectrum of the emitted light (color)
- Stability of intensity
- Handling
- Lifetime

For offering the optimal solutions for all these applications and requirements we put together the following selection of standard light sources:

LED light sources:

30-70 W with various color temperatures

Halogen fiber light sources:

30 W, 100 W and 150 W, with continuous light regulation, color temperatures of 3,000-3,400 K (white light)

Metal halide fiber light sources:

100 W and 150 W, with continuous light regulation, color temperatures of 4,200 and 5,200 K

To guarantee the optimum light transfer to the object, we offer two types of light guides:

Light guides with glass fiber bundles:

These consist of many highly flexible, individual glass fibers that are optically processed, suitable for low and medium outputs in the visible spectrum. Their transmission is approx. 50 %. The following versions are available:

The following options are available:

- Flexible, protected with PVC or metal tubing
- Flex and stay metal coil (swan neck)
- One, two, three or more arms
- Ring-shaped or inline
- Customer-specified light guide

Liquid fiber light guides:

Flexible, liquid-filled PVC tubing that is sealed on the ends with quartz windows. This type of fiber light guide is ideal especially for applications in which high transmissions (approx. 85 %) and high powers are required. The three standard liquid fiber light guides cover a wavelength range from 300 nm (UV) to 2,000 nm (IR).

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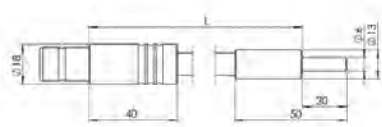
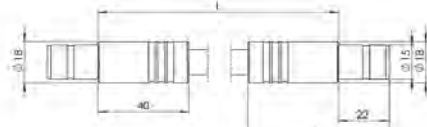
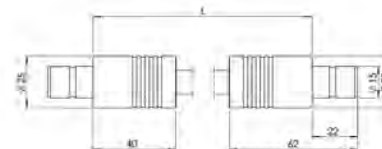
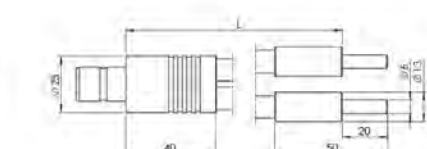
Dimensions Flexible Light Guides

Flexible Fiber Light Guide

Flexible light guides are an easy-to-use and flexible source of pinpoint illumination.

Qioptiq offers flexible glass fiber light guides in metal tubing as standard products in single, dual and triple arm models. These light guides are equipped with the Light Source Connector LA and metal ferrules (black-anodized aluminum).

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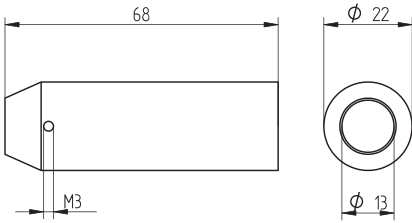
Focusing Lens

- For magnifying the operating distance
- For positioning on the end of the light guide of the series SL, SLZ and FL
- Fixed in place by a locking screw

- Weight: 0,032 kg
- Diameter: 22 mm
- Length: 68 mm
- Angle of Aperture (2Φ/2): bis zu 20°
- Clamp diameter (Fiber Ferrule): 13 mm
- Material: Black anodized aluminum and glass

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Produktbezeichnung	Artikel-Nr.
<i>Item Title</i>	<i>Part No.</i>
Focusing lens	G405815000

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Fiber optic ring light with an inner diameter of 66 mm for usage with a cold light source.

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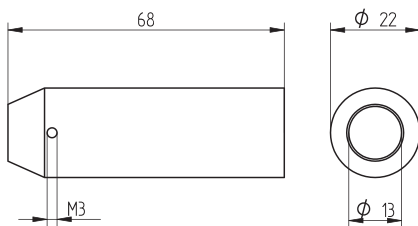
Focusing lens

- For magnifying the operating distance
- For positioning on the end of the light guide
- Fixed in place by a locking screw
- Weight: 0.032 kg
- Diameter: 22 mm
- Length: 68 mm
- Angle of Aperture (2Φ/2): bis zu 20°
- Clamp diameter (Fiber Ferrule): 13 mm
- Material: Black anodized aluminum and glass

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Produktbezeichnung	Artikel-Nr.
Item Title	Part No.
Focusing lens	G405656000

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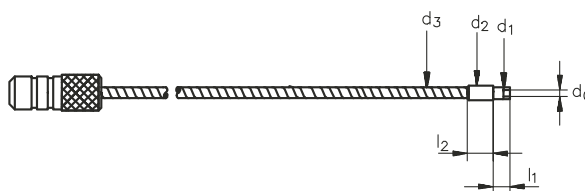
Liquid Fiber Light Guides LLG VIS

Single arm, liquid-filled light guide jacketed with black PVC with light guide connection LA. The light exit is equipped with an aluminium ferrule. The ends of the light guide are sealed through fused silica windows.

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- Can be used with all Qioptiq fiber light sources
- Numerical aperture (2α) of 72° is considerably higher than the conventional glass fiber light guide.
- Wavelength Range: 340-800 nm
- Max. Transmission (380-720 nm) > 70 %
- Temperature range - 5 °C to + 35 °C

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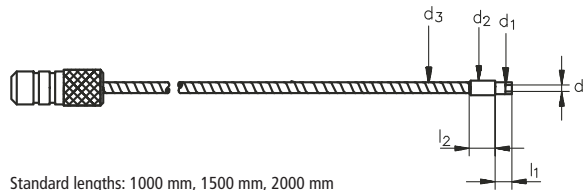
Liquid Fiber Light Guides LLG IR

Broadband, single armed light guide like LLG VIS, but optimized for the VIS and IR wavelength range.

- Can be used with all Qioptiq fiber light sources
- Numerical aperture (2α) of 72° is considerably higher than the conventional glass fiber light guide.
- Wavelength Range: 420-2000 nm
- Max. Transmission (420-1400 nm) > 70 %
- Temperature range - 5 °C to + 35 °C

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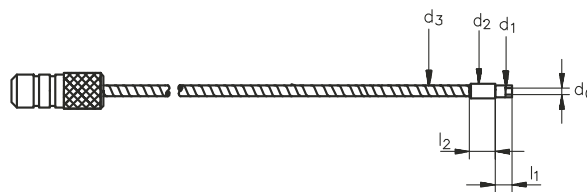
Single arm light guide like LLG VIS for the transmission range of 300 to 650 nm.

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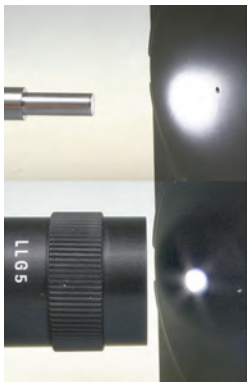
Focusing Lens Attachments FA-LLG

Light emission of a Liquid Fiber Light Guide without (top) and with focusing adapter (bottom)

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Focusing Lens Attachments FA-LLG

Attachment for all liquid fiber light guides with 3 and 5 mm diameters.

Focuses the divergent light emitted by the fiber light guide to a point that has a diameter of approx. 5 mm.

The 3-lens quartz glass system allows the attachment to be used in the spectral region from UV to near IR. The focusing lens attachment mounts directly to the end of the liquid fiber light guide LLG.

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Cold Light Sources CLS-LED3

LED Cold Light Sources by StarLight are of high quality. Their sophisticated design and the high-class LED ensure reliability, efficiency and durability.

The Cold Light Source LED3 is passively cooled (no fan) which makes it perfect for clean room applications. Its luminous intensity exceeds a conventional 150 W Halogen Cold Light Source.

There are various versions of this fan-less cooled LED Cold Light Source. You can choose between high lumen output and high color rendering index. Also there are models with natural white and warm white light. So you can find the right LED Cold Light Source for your application.


- Flicker-free light by optimized electronic
- Robust aluminum body with modern design
- Adaptable for several light guides (adapters optionally available)
- Horizontal vertical positioning, stackable
- Illuminant: High-Power-LED 30 W
- Input voltage of cold light source: 15 V DC
- Max. power consumption: ca. 36 W
- Ø light guide connection: 15 mm (changeable with adapter)
- LED life span: 50.000 h (manufacturer specification)
- Material: anodized aluminum
- Color: titanium, grey
- Weight: ca. 3,7 kg
- Input power supply: (EU|UK|US) 100-240 V AC, 50/60 Hz
- Output power supply: 15 V DC, 2.400 mA
- Dimensions (L×W×H): ca. 260 × 175 × 80 mm

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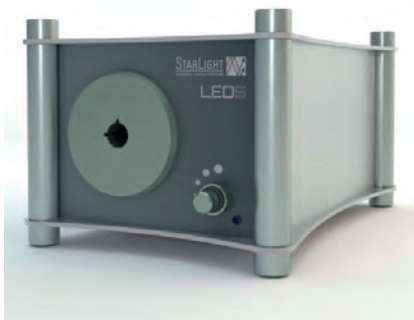Cold Light Source CLS-LED5


LED Cold Light Sources by StarLight are of high quality. Their sophisticated design and the high-class LED ensure reliability, efficiency and durability.

The actively cooled LED Cold Light Source LED5 is a true all-rounder with an excellent price-performance ratio. The great advantage of this Cold Light Source is the high light yield for light guides with a small diameter as well as for those with a large diameter.

- Flicker-free light by optimized electronic
 - Glare protection by light guide detection
 - Robust aluminum housing with modern design
 - Adaptable to all kind of light guides (adapters optionally available)
 - Horizontal and vertical positioning, stackable
 - Build-in power supply
 - Power cable included
- Illuminant: High-Power-LED 50 W
 - Luminous flux with light guide 1 m × Ø 5 mm: 400 lm
 - Luminous flux with light guide 1 m × Ø 9 mm: 810 lm
 - Luminous flux with light guide 1 m × Ø 13 mm: 980 lm
 - Color temperature: 7.700 K
 - Color rendering index: 70
 - Max. power consumption: 62 W
 - Operating voltage: 100-240 V AC, 50/60 Hz
 - Ø light guide connection: 15 mm (changeable with adapter)
 - Active diameter light guide: max. 13 mm
 - Ambient temperature: 5°C to 35°C
 - LED life span: 50.000 h (manufacturer specification)
 - Width: 174 mm
 - Height: 137 mm
 - Length: 259 mm
 - Material: anodized aluminum
 - Color: titanium, grey
 - Weight: ca. 2,7 kg
 - Diameter light guide connector: 15 mm
 - Light color: cold white
 - Max. power consumption: 62 W
 - Supply voltage: 100-240 V AC, 50/60 Hz
 - Dimensions (L×W×H): ca. 260 × 175 × 140 mm

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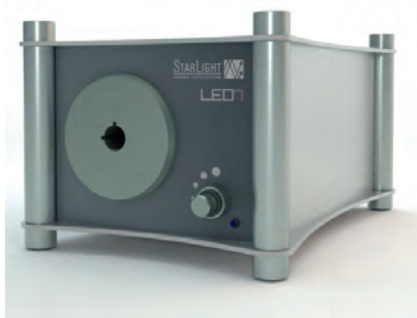Cold Light Sources CLS-LED7

LED Cold Light Sources by StarLight are of high quality. Their sophisticated design and the high-class LED ensure reliability, efficiency and durability.

The actively cooled LED Cold Light Source LED7 is an uncompromising power pack. It is the perfect choice for applications where extreme brightness is needed. The low-noise fan provides a long lifespan of the LED, even in permanent use.

- Flicker-free light by optimized electronic
- Glare protection by light guide detection
- Robust aluminum housing with modern design
- Adaptable to all kind of light guides (adapters optionally available)
- Horizontal and vertical positioning, stackable
- Build-in power supply
- Power cable included
- Illuminant: High-Power-LED 70 W
- Luminous flux with light guide 1 m × Ø 9 mm: 1350 lm
- Color temperature: 6.900 K
- Color rendering index: 75
- Max. power consumption: 78 W
- Operating voltage: 100 - 240 V AC, 50/60 Hz
- Ø light guide connection: 15 mm (changeable with adapter)
- Ambient temperature: 5°C to 35°C
- LED life span: 50.000 h (manufacturer specification)
- Width: 174 mm
- Height: 137 mm
- Length: 259 mm
- Material: anodized aluminum
- Color: titanium, grey
- Weight: ca. 2,7 kg
- Diameter light guide connector: 15 mm
- Light: color cold white
- Supply voltage: 100-240 V AC, 50/60 Hz
- Dimensions (L×W×H): ca. 260 × 175 × 140 mm

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IL100-c lamp

Mounting modules to IL100-c

LED Industrieleuchte IL100-c LED-Industry Lamp IL100-c

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LED-Industry Lamp IL100-c

The LED Machine Lamp IL100-c is a versatile and robust LED Lamp for industrial applications with changing requirements. Typical scope of this Machine Lamp you can find in printing, packaging, tool manufacturing, metalworking, automation, machine vision and measuring.

The eco-friendly and durable LED technology in combination with the optimized electronic enables a safe and easy usage.

Thanks to the innovative concept "modul-LED" you can change the LED module fast and easy, even during operation. So you can choose between two light colors.

With optionally available accessories you can mount the LED Machine Lamp IL100-c wherever you want and adjust the light spot on the area to be illuminated.

The robust anodized aluminum body and the protective glass avoid mechanical or chemical damage of the LED module. Furthermore, dust or water can't come into the lamp. So the LED Machine Lamp is suitable for applications in harsh environments.

- *Insensitive to vibrations*
- *Fast and simple replacement of the LED module*
- *No generation of heat in light direction*
- *Homogenous and flicker-free illumination*
- *Type of LED: 3 W Power-LED*
- *Illuminant: LED module 21 mm (exclusive)*
- *Connection cable with industrial connector (5-pol)*
- *Dimensions table control (LxWxH): 78 mm x 38 mm x 21 mm*
- *Length of feed line: 3 m*
- *Ø LED module: 21 mm*
- *Diameter: 25 mm*
- *Length: 96,9 mm (head with protection cap)*
- *Material: anodized aluminum*
- *Color: titanium*
- *Input power supply: (EU|UK|US) 100-240 V DC, 50/60 Hz output power supply 24 V DC, 1 A*
- *Cable length power supply: ca. 1,4 m*
- *Dimensions with package: 255 mm x 155 mm x 90 mm*
- *Weight with package: approx. 540 g (with power supply, TC and LED module)*
- *Weight controller: 50 g*

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Modules for IL100-c

The innovative concept "modul-LED" allows a fast exchange of LED modules. There are two versions of the LED-module differing in light color.

Modules for IL100-c

Module für IL100-c **Modules for IL100-c**

Produktbezeichnung	Farbtemperatur (K)	Artikel-Nr.
<i>Item Title</i>	<i>Color Temperature (K)</i>	<i>Part No.</i>
Module for IL100-c, Flood 25°-6000K	6000	G405818000
Module for IL100-c, Flood 25°-4000K	4000	G405819000

NEW

Mikrobankbeleuchtung LED ML

Microbench and Microscope light source ML3

Mounting example of ML3 to Microbenchsystem

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Microbench Light Source LED ML

The super-bright Microbench lights in the LED ML Series emit homogenous and flicker-free light. A digital potentiometer is used for processor-controlled adjustment of the brightness with no noticeable change in the color temperature. The most recent brightness setting can be automatically stored in memory for the next use.

The LED ML Series Microbench Lights offer you all the advantages of the latest in LED technology, as an energy efficient and environmentally friendly alternative to halogen lights. The high-quality 3 W power LED has a service life of approximately 50,000 hours. This ensures almost six years of continuous operation without complicated, expensive bulb replacements.

- Homogeneous and flicker-free illumination
- Energy-saving, environmentally friendly lighting technique
- Color LEDs in red, blue and green
- White light in pure-white, natural-white and warm-white
- UV-LED with 405 nm, 375 nm and 365 nm
- Digital brightness control; no noticeable change in color temperature
- Most recent brightness setting stored in memory
- Long LED Service Life: Approx. 50,000 h
- Light: 3 W power LED
- LED Service Life: Approx. 50,000 h
- Weight: 0.1 kg
- Dimensions: Length: 120 mm (4.72 in.), Ø: 30 mm (1.18 in.)
- Housing Material: Aluminum, black anodized
- Operating Voltage: 12 V DC
- Operating Voltage, Power Supply: 100-240 V AC, 50-60 Hz
- Equipment Included: 24 V DC power supply

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X-Cite® Fluorescence Illumination Solutions

X-cite products are originally developed by Lumen Dynamics, a leader and volume manufacturer of lamp and LED-based UV curing and fluorescence illumination systems. Since 2013 Lumen Dynamics is part of Excelitas Technologies.

More details of our famous x-Cite illumination products are shown under www.ldgi.com/de/x-cite/

X-Cite® Product Overview

X-Cite® offers the Life Science and Analytical Instrumentation market a broad range of innovative fluorescence illumination and measurement solutions with unsurpassed reliability and control capabilities. Whether observing fixed or live cells, X-Cite® offers a complete range of lamp and LED microscope light sources that optimize imaging and ensure greater data reliability. These X-Cite® products are known for maximum stability and superior illumination uniformity for fluorescence microscopy applications, with lamp-based solutions that provide the convenience of pre-aligned, long-life lamps and easy installation. Delivering high intensity LED illumination, the X-Cite® XLED1, X-Cite® 120LEDBOOST, X-Cite® 110LED, and X-Cite® TURBO feature intuitive controls and no bulbs to install, making set-up and operation simpler than ever.

Fluorescence Microscope Light Sources

X-Cite® offers a wide range of lamp and LED light sources and fluorescence illumination systems ideal for a variety of microscopy applications. From our standard X-Cite® 120Q model used for routine assays, to our most advanced high-end X-Cite® XLED1 and X-Cite® TURBO, X-Cite® has the right fluorescence illumination system for your application.

Power Measurement in Fluorescence Imaging

The X-Cite® Power Measurement System measures optical power where it matters most — at the objective plane. With consistent illumination levels, reproducible imaging results can be ensured while increasing the integrity of your research.

X-Cite® Lamp Systems Overview

Fluorescence Illumination - In Control

	X-Cite 120Q	X-Cite 120PCQ	X-Cite 200DC	X-Cite exacte
Lamp/LED Type	120 W AC Lamp	120 W AC Lamp	200 W DC Lamp	200 W DC Lamp
Wavelength Range	320-750 nm	320-750 nm	340-800 nm	340-675 nm
Lamp/LED Warranty	2,000 hours	2,000 hours	2,000 hours	2,000 hours
Mercury Content	20 mg	20 mg	30 mg	30 mg
Intensity Control	5-step (optional)	5-step	0-100%	100-step
Shutter Speed	N/A	~200 ms	5-6 ms (via TTL)	5-6 ms (via TTL)
PC Control	N/A	Yes	TTL shutter	Yes
LLG Detection	No	No	Yes	Yes
Consumables	Bulb, liquid light guide	Bulb, liquid light guide	Bulb, liquid light guide	Bulb, liquid light guide
Coupling	Light guide	Light guide	Light guide	Light guide
Imaging Requirements	<ul style="list-style-type: none"> High illumination power Medium term stability¹ 	<ul style="list-style-type: none"> PC control Automated imaging 	<ul style="list-style-type: none"> Short exposure times² High speed shuttering 	<ul style="list-style-type: none"> PC control Automated imaging Multi-day experiments

¹Medium term = exposures greater than 100 ms

²Short term = exposures less than 50 ms

X-Cite® LED Systems Overview

Fluorescence Illumination - In Control

	X-Cite 110LED	X-Cite 120LEDBoost	X-Cite 120LEDmini	X-Cite XLED1	X-Cite Turbo
LED Type	Broadband	Broadband	Broadband	4 switchable wavelengths	6 switchable wavelengths
Wavelength Range	360-660 nm	370-700 nm	370-700 nm	10 LED options (360-750 nm)	380-650 nm
LED Warranty	25,000 hours	25,000 hours	20,000 hours	20,000 hours	15,000 hours
Mercury Content	None	None	None	None	None
Intensity Control	1% increments	1% increments	1% increments	0.1% increments	1% increments
Analog Intensity Control	No	No	No	No	Yes
Shutter Speed	1 ms USB 100 µs TTL	500 µs USB 50 µs TTL	1 ms USB 85 µs TTL	500 µs USB 10 µs TTL	1ms USB 100 µs TTL
PC Control	Yes	Yes	Yes	Yes	Yes
Consumables	Liquid light guide	None	None	Liquid light guide	Liquid light guide
Coupling	Light guide	Direct coupling	Direct coupling	Light guide	Light guide
Imaging Requirements		<ul style="list-style-type: none"> Broadband light Stability Minimal sample exposure 		<ul style="list-style-type: none"> Fast wavelength switching Pulsing to increase cell viability Stable and repeatable Advanced automation 	<ul style="list-style-type: none"> Turbo power in 500-600 nm region Fast wavelength switching Stable and repeatable

X-CITE® Power Measurement and Accessories

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LED coupler

- *To increase coupling efficiency into glass fiber bundles*
- *Length: 12.30 mm (optics), 9.70 mm (mechanics)*
- *Diameter: M12x0.5*
- *Wavelength: 400 – 650 nm*
- *Transmission: > 96 %*
- *Captured beam angle of LED: $\pm 45^\circ$*
- *Coupling efficiency: > 60%*
- *LED diameter: 3 mm*
- *Distance to LED X: 0.50 mm*
- *Fiber diameter: 5 mm*
- *Distance to fiber Y: 3.50 mm*
- *Temperature range: 0°- 80°C*

Index

[Alphabetical and Numerical Index](#)
[Terms and Conditions](#)

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616
628
640

Content

Alphabetical Index
Numerical Index
Terms and Conditions

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Alphabetical Index

A

Achromatic Retarders	402
Achromats VIS, Positive; Dia. 3 mm to 31.5 mm, Mounted	290
Achromats VIS, Positive; Dia. 3 mm to 31.5 mm, Unmounted	288
Achromats VIS, Positive; From 40 mm Diameter, Unmounted	292
Adapter 1" to W0.8" or Ø 25 mm	62, 122
Adapter 1.035"-40UN / 30	63
Adapter 16	64
Adapter 16/25	58
Adapter 19	64
Adapter 23.2	65
Adapter 23.2 to Ø 25	66
Adapter 25/15	66
Adapter 25/C-Mount	61
Adapter 30/25	59
Adapter 30/C-Mount	61
Adapter 35/30	60
Adapter 35/30-3.5°	60
Adapter Bench Systems 30 & 1.035"-40 UN	63
Adapter C-Mount/Microbench	62
Adapter Collar	59
Adapter FIC / Microbench	576
Adapter M35x0.5/30	60
Adapter Plate	68
Adapter Plate MF-25	235
Adapter Ring C	123
Adapter Ring C-W0.8	123
Adapter W0.8" to 1"	65, 122
Adapter W0.8"x1/36	65
adjust.x Medium L-Type	234
adjust.x Small L-Type	233
Adjustable Height Column 75	200
Adjustable Holder 25	32
Adjustable Holder 35-M	194
Adjustable Holder 55-M	194
Adjustable Mirror Insert 30	51
Adjustable Optic Holder 31.5	56
Adjustable Prism Insert 30	50
Adjusting Tube C30 with Locking Collar	118
Adjustment Plate N	94
Allen Wrench Set	213
Aluminized Plane Mirrors - Circular plane	342
Aluminized Plane Mirrors - Elliptical plane	344
Aluminized Plane Mirrors - Rectangular plane	341
Amplifier for Phase Modulators	541
Angle Bracket 40	204
Angle Bracket 80-M	205
Aperture Set, Filter Mount and Adjustable Iris Diaphragm	467
Apo-Rodagon D	438
Apo-Rodagon HR	435
Apo-Rodagon-N	439
Aspheric Condenser Lenses, Crown-glass	275
A-Zoom Micro	472
A-Zoom2 10X Single-Objective Microscope	470
A-Zoom2 40X Single-Objective Microscope	471

B

Base Plate 80x40-M	207
Base Plate N	91
Base Plate N 40	92
Base Plate N 50	92
BBO Double Pockels Cells DBBPC Series	553
BBO Pockels Cells BBPC Series	552
BBO Pockels Cells for High Switching Rates DBBPC HR Series	554
Beam Expanders, Variable Magnification	416

Beam Blocking Shutter, Mechanical	576
Beam Deflector Cube C	109
Beam Deflector Cube C-5	110
Beam Dump	116
Beam Expanders, Motorized	418
Beam Steering Mirror Assembly	54
Beam-Deflector Cube	36
Beamsplitter Cube Insert, Adjustable	112
Beamsplitter Cubes NIR	322
Beamsplitter Cubes UV	320
Beamsplitter Cubes VIS	321
Beamsplitter Plates IR	327
Beamsplitter Plates NIR	326
Beamsplitter Plates UV	323
Beamsplitter Plates VIS-NIR	324
Beamsplitter Plates with Different Splitting Ratios	328
Beamsteering Mirror Mount, adjustable	111
Blue Colored Glass Filters	309
Broadband Anti-Reflective Coating ARB 1	369
Broadband Anti-Reflective Coating ARB 1 IR	369
Broadband Anti-Reflective Coating ARB 1 UV	369
Broadband Anti-Reflective Coating ARB 2 IR	371
Broadband Anti-Reflective Coating ARB 2 NIR	370
Broadband Anti-Reflective Coating ARB 2 UV	370
Broadband Anti-Reflective Coating ARB 2 VIS	370
Broadband Anti-Reflective Coating ARB 3	371
Broadband Anti-Reflective Coating ARB 3 NIR	371
Broadband Anti-Reflective Coating ARB 4	372
Broadband Anti-Reflective Coating ARB 5	372
Broadband Beamsplitter Coating TBP VIS	384
Broadband Beamsplitter Coating TBP VIS NIR	384
Broadband Hybrid Beam-splitter Coating TBW-H VIS	386
Broadband Laser Mirror Coating DLB VIS	382
Broadband Mirror Coating DLB UV	381
Broadband Mirror Coating DLB SBB	382
Broadband Narrowband Anti-Reflective Coating ARBS	373

C

Cap 30/25	59
Carrier 65	175
Carrier FLR 40	172
Carrier Positioner X 95 - 30	181
Carrier X 95 - MB	182
Carriers FLR 95	178
Carriers X 95	180
Centering Lens Holder 100	198
Centering Mounting Plate/Holder 25	40
CF 30 Controller for Piezo Electric Inertial Motors	161
Circlip Kit	75
Clamp Holder for Prism	55
Clamping Fork	199
Clamping Set for Stages	151
Clamps SP 4	212
Clamps SP 4 and SP 4-10	183
Closing Caps	72
Closing Tube C30x6	107
C-Mount Adapter for HeNe Lasers with Outer Diameter of 45 mm	574
C-Mount Adapter for HeNe-Laser with Outer Diameter of 35 mm	574
C-Mount Adapter M17	121
C-Mount Adapters for Compact Objectives	448
C-Mount Microbench Adapter	123
C-Mount-Adapter-M17	61
CN 30 Controller for Piezo Electric Inertial Motors	161
Cold Light Mirror KS 93 / 45°	318
Cold Light Source CLS-LED5	605
Cold Light Sources CLS-LED3	604
Cold Light Sources CLS-LED7	606

Collar N	100
Collimator Holder	71
Color Filters DC	314
Color Filters DT	312
Colored Glass Filters, Different Edge-Positions	311
Compact Objectives	444
Compression Spring, 60x8x0.7	50
Condenser C35 with Ancillary Lens	584
Condenser Combination, $f=16$ mm, Crown Glass	293
Condenser; One-Lens 220 mm to ∞ ; UV	596
Condenser; One-Lens 220 mm to ∞ ; VIS	596
Condenser; Two-Lens; 105-385 mm; UV	595
Condenser; Two-Lens; 105-385 mm; VIS	595
Connecting Clamp, 14 mm Diameter	201
Connecting Tube C30 FC / C30 FSMA	119
Connecting Tube C30 with Adjustable Mount 6	120
Connection cables for piezo positioners	161
Connection Clamp 14 V	202
Connector 80 V	37
Corner Connector 40/80/40 with Threads	36
Corner Connector T	37
Counter Support	49
Cross Carrier X 95	182
Cube	35
Cube N	93
Customized Micro Objectives	481
Cuvette Holder	71
Cylindrical HeNe Lasers w/o Power Supply, Outer Diameter 31.7 and 44.25 mm	572
Cylindrical HeNe Lasers w/o Power Supply, with Shutter, Outer Diameter 35 and 45 mm	571

D

Dielectric Broadband Beamsplitter Coating TBW UV	385
Dielectric Broadband Beamsplitter Coating TBW VIS	385
Dielectric Coated Plane Mirror DLHS 755-840 nm, Ti:Sa	351
Dielectric-Coated Plane Mirrors DLB SBB 350-950 nm	354
Dielectric-Coated Plane Mirrors DLB UV 230-400 nm	353
Dielectric-Coated Plane Mirrors DLHS 532 nm	349
Dielectric-Coated Plane Mirrors DLHS 633 nm	350
Dielectric-Coated Plane Mirrors DLHS NIR 1064 nm	352
Dielectric-Coated Plane Mirrors DLHS NIR 1540 nm	352
Dielectric-Coated Plane Mirrors DLHS UV 248 nm	346
Dielectric-Coated Plane Mirrors DLHS UV 266 nm	346
Dielectric-Coated Plane Mirrors DLHS UV 308 nm	347
Dielectric-Coated Plane Mirrors DLHS UV 351-355 nm	348
Dielectric-Coated Plane Mirrors DLHS UV 405 nm	349
Dielectric-Coated Plane Mirrors DLHS VUV 157 nm	345
Dielectric-Coated Plane Mirrors LBSM-NIR	356
Dielectric-Coated Plane Mirrors LBSM-UV	355
Dielectric-Coated Plane Mirrors LBSM-VIS	355
Dielectric-Coated Plane Mirrors LBSM-YAG	356
Digital Pulse Amplifier DIV 20	540
Diode and DPSS Laser	562
Dispersion Plate in Retaining Mount	584
Double Mirror Coating DLHD UV/VIS	380
Double Mirror Coating DLHD VIS/NIR	381

E

Electrical Lens Driver	281
Electrically Focus Variable Lens EL-10-30	278
Electrically Focus Variable Lens EL-16-40	280
Eyepiece Adapter 23.2/30	66
Eyepiece Adapter C30	120

F

Faraday Isolators, Overview	511
Faraday Isolators, Single Stage	512
Faraday Isolators, Two Stage	525
FC Fiber Adapter 3.5°	67
FC/FSMA Fiber Adapter	67
Fetura+	464
Fiber Illumination System	597
Fiber Illumination Systems, Overview	597
Fiber Light Guide, Flexible	599
Fiber Optic Adapter N FC	98
Fiber Optic Adapter N FSMA	99
Fiber Optic Ring Light	599
Filter Support C	114
Fine Adjustment C30	119
Fine-Adjustment Screws, Flat Bushing	159
FishEye Objective 190°	445
Flat Rails 65	174
Flat Wrench	125
fle.X-Plate	42
Flexible Fiber Light Guide	598
FLS 95 Rail System	179
Fluorescence Microscope Light Sources	610
Fluorescent Cube Holder	467
Focusing Collimators	73
Focusing Lens	598
Focusing lens	600
Focusing Lens Attachments FA-LLG	603
Focusing Mount C30	118
Focusing Mount C35	120
Focus-Variable Lenses from Optotune	276
Frame Holder 100-M	199
Front Surface Mirror Silver Coating RAGV	377
Front Surface Mirrors for Illumination Systems	358
F-Theta-Ronar Lenses	408
Fused Silica Brewster Plates	303
Fused Silica Brewster Polarizer	394

G

Gimbal Mirror Holder MA	236
Glan-Taylor Prism Polarizers	392
Glan-Thompson Prism Polarizers	393
Glass Manager	496
Gold Plane Mirrors, Laser Quality	334
Goniometer Stages 60-60 / 60-80	152
Green Colored Glass Filters	310

H

Halogen Lamp Power Supply PSH 100	590
Hard Seats	232
HD Micro Objectives	294
Heat-Absorbing Filters	316
Heat-Protection Filters CALFLEX™ 3000 SP	317
Heat-Protection Filters CALFLEX™ B1/K1	317
Heat-Protection Filters CALFLEX™ C	318
HeNe Laser and Accessories	568
HeNe Lasers, Power Supply	573
High-Power Doubleband Anti-Reflective Coating ARHDS UV-VIS	374
High-Power Doubleband Anti-Reflective Coating ARHDS VIS-NIR	374
High-power laser mirror coating DLHS Ti:Sa	380
High-Power Laser Mirror Coating DLHS VIS	379
High-power Narrowband Anti-Reflective Coating ARHS	373

Alphabetical Index

High-Power Narrowband Anti-Reflective Coating ARHS UV	373
Holder 150	221
Holder 25/30/35/M35x0.5	32
Holder for Micrometer	49
Holder N 20	90
Holder N 40	91

I

iFLEX-Agile Series cw-OPO	577
iFLEX-iRIS Accessories	567
iFLEX-iRIS CDRH Compliant Interlock	566
iFLEX-iRIS Fiber Coupled Laser	564
iFLEX-iRIS Laser	562
iFLEX-iRIS OEM power supply	567
Incident Light Microscope Objectives	293
Infinity Macro Lens or Wide Field Lens	468
Inspec.x L 4/105	432
Inspec.x L 4/60 and 4/100	434
Inspec.x L 5.6/105	430
Inspec.x L 5.6/105 float	431
Inspec.x L Prism Module	433
Inspec.x M	429
Instrument Flange	68
IR Broadband Beamsplitter Coating TBP IR	385
IR High-Power Laser Mirror Coating DLHS NIR	380
Iris Diaphragm 50 in Mounting Ring	197
Iris Diaphragm Nanobench	214
Iris Diaphragm Set	116
Iris Diaphragm Tube Mounting System C	215
Iris Diaphragms Microbench	214
Iris Diaphragms with Different Outer Diameters	216
Iris Diaphragms, Completely Closing	216
Isolator with 8 mm Aperture, SI Series	518
Isolators with 2 mm Aperture, SV Series	513
Isolators with 3.5 and 5 mm Aperture, SC Series	514
Isolators with 5 mm Aperture, SV Series	517
Isolators with a Broad Tuning Range	521
Isolators with Magneto-Optical Crystal Film	520
Isolators, Tunable	522

K

KD*P Brewster Pockels Cell BPC 8	552
KD*P Double Pockels Cells DPZ Series	551
KD*P Pockels Cells CIQS Series	549
KD*P Pockels Cells CPC Series	548
KD*P Pockels Cells LM Series	547
KD*P Pockels Cells SPC 4 Series	550
Knurled Ring C30x8	122

L

Lamphouse IL	586
Laser Adapter C 35 for HeNe Laser with Outer Diameter 31.7 mm	575
Laser Beam Expander System 10x	420
Laser Beam Expander System 4x and 7x	420
Laser Modulators LM 0202	538
Laser Modulators LM 13	536
Laser Speckle Reducer	422
Laserset	25
LED Coupler	612
LED Light Sources	604
LED Light Sources HighLED	583
LED-Industry Lamp IL100-c	607
Lens Design Elements Poster	502
Lens Mount	222

Lens mounts for diameter 3 mm - 31.5 mm	219
Lens Protection Cap	222
Lens Tube C35x50	108
Lens Turret 4x with C-Mount Adapter	74
Lid C	110
Light Trap	116
Lightweight Breadboards CleanTop® 75 Series	208
Linear Positioners TM / TMS	134
Linear Stages for Microbench	45
Linear Stages T 10 M/F and TS 10 M/F	140
Linear Stages TB	136
Linear Stages TL 45 / TL 65	142
Linear Stages TL 45 / TL 65 with Digital Micrometer	143
Linear Stages TL 65 with Side Drive	144
LINOS Focus-Ronar Lenses 1064 nm/532 nm	415
LINOS Focus-Ronar Lenses, 355 nm, Fused Silica	414
LINOS F-Theta-Ronar Lenses 1064/1030-1080 nm	411
LINOS F-Theta-Ronar Lenses 1550/1900-2000 nm	412
LINOS F-Theta-Ronar Lenses 340-360 nm	408
LINOS F-Theta-Ronar Lenses 515-540/532 nm	409
LINOS F-Theta-Ronar Lenses 940-980 nm	410
LINOS Motorized Beam Expanders	418
LINOS Variable Magnification Beam Expanders	416
Liquid Fiber Light Guides LLG IR	601
Liquid Fiber Light Guides LLG UV	602
Liquid Fiber Light Guides LLG VIS	600
LM1 Optic Mount	226
LM2 Optic Mount, Compact	228
Low Outgassing Isolators with 3.5 mm aperture, LO Series	516
LWD Objectives	463

M

MachVis Software	500
MachVis Workflow	501
Macro CCD Lenses	442
mag.x System 125	452
Magnetic Base	207
Magnetic Cage Assembly Swapper Plates	31
Magnetic Clamp 50-M6 (4 ea.)	208
Material Editor	498
Mechanical Accessories	446
Mercury Arc Lamp	588
Mercury Arc Lamp Power Supply PSQ 50AC	592
MeVis-C	426
MeVis-C/CF Traffic	428
MeVis-CF	427
Micro Optical Customized Components	481
Micro Optics, General	480
Microbench	11
Microbench - Nanobench Adapter	58
Microbench Adapter	98
Microbench Adapter for HeNe-Laser with 31.7 mm Outer Diameter	575
Microbench Adapter IL	594
Microbench Alignment Plate	75
Microbench Basic Set	22
Microbench Light Source LED ML	609
Microbench Light Sources	582
Microbench Mechanics Test Kit	26
Microbench Set Mechanics	19
Microbench Set Optics	21
Microbench Set O-UV	24
Micrometers G	158
Mirror Holder 45°	95
Mirror Holder 45°/7x12	96
Mirror Mount MF-25PF	235
Modules for IL100-c	608
Motorized Rotation Mount	48
Mount 100	221

Mount 31.5	117
Mounted Pinholes	217
Mounting Base	206
Mounting Base 80x15	207
Mounting Bases	34
Mounting Bracket 40	39
Mounting Bracket 40/65	594
Mounting Bracket WM 50-1/80-1 for Linear Stages TB	138
Mounting Bracket WM 50-2/80-2 for Linear Stages TB	139
Mounting Column 14	200
Mounting Components	145
Mounting Plate 220x140-M	210
Mounting Plate 220x40 mm	210
Mounting Plate 25 - T10	30
Mounting Plate 25 with Mounting Holes	28
Mounting Plate 25/30, Latchable, 10°	29
Mounting Plate 25/30/35, Latchable	28
Mounting Plate 25E/30E	30
Mounting Plate 80x115-M	211
Mounting Plate 80x80	210
Mounting Plate FLS	72
Mounting Plate M35x0.5, also Latchable	31
Mounting Plate MP 80.U	211
Mounting Plate N	89
Mounting Plate N, latchable and rotatable	90
Mounting Plates	209
Mounting plates 25	209
Mounting Plates 25/30/35/1"	27
Mounting Ring for Microbench	196
Mounting Rings for Mounted Optics	196
Mounting Tool for Beamsplitter Cubes	52
Mounts for Post Use	202
MS 30 Miniature Translation Stage	162
MS 38 Miniature Translation Stages with Free Opening	163
MT 25 Kinematic Mirror Tilting Stage	165
MX 25/35 Miniature XYZ Positioners with Piezo Electric Inertial Drive	164

N

Nanobench	79
Nanobench Basic Set	88
Nanobench Mechanics Test Kit	84
Nanobench Set Mechanics	85
Nanobench Set Optics	87
N-BK7 Plane Plates	300
Neutral Wedges	316
Neutral-Density Filters	314
NIR Broadband Beamsplitter Coating TBP NIR	384
NIR-Blocking Filters CALFLEX™ X	319

O

Object Holder 50-M	198
Optem Camera Mounts	459
Optem Camera Tubes	460
Optem Fusion Basic Modules	459
Optem Fusion Illumination	468
Optem Fusion Motorization	469
Optem Lower Function Modules	462
Optem Lower Lens	462
Optem Zoom Modules	461
Optic Holder 63-M	195
Optic Mount 1", C30x15	117
Optic Mount 63 (40, 50, 50.8 mm)	220
Optic Support 25	54
Optical Contacted Polarizing Beamsplitter Cubes	398
Overview Table Adapter	57

P

Partially Reflective Laser Mirror Coating TDLS	383
Penta Prisms	306
Phase Modulator PM 25	534
Phase Modulator PM-C-BB	535
Pinholes for Spatial Filters	218
Plane Mirror SILFLEX	357
Plano-Concave Lenses, Mounted (Fused Silica)	270
Plano-Concave Lenses, Mounted (N-BK7)	269
Plano-Concave Lenses, Unmounted (Fused Silica)	268
Plano-Concave Lenses, Unmounted (N-BK7)	267
Plano-Convex Lenses, Mounted	259
Plano-Convex Lenses, Mounted (Fused Silica)	260
Plano-Convex Lenses, Unmounted	256
Plano-Convex Lenses, Unmounted (Fused Silica)	258
Plate Holder	197
Pockels Cells Positioner	556
Pockels Cells Product Overview	546
Polarization Filter 80 in Rotary Mount	404
Polarization Filter in Retaining Mount	585
Polarization Filter Type VIS 4 K	403
Polarizing Beamsplitter Cube Coating TSWP	386
Polarizing Beamsplitter Cubes	395
Polyamide Plugs	78
Porro Prisms, 90° (Isosceles)	306
Positioning Ring 30	53
Post 25-200 with Clamp	203
Post 38	204
Power Measurement in Fluorescence Imaging	610
Power Supply for Cylindrical HeNe Lasers	573
Precision Adjustment Screws G, bushing with connection thread	158
Precision Lead Screws Lees	160
PreDesigner	499
Prism Holder	96
Prism Stages 40 S / 60 S	157
Prism Support 25	55
Prism Table 25	56
Prism Table 25, Adjustable	52
Prism Table Insert, Adjustable	111
Prisms, 60° (Equilateral)	307
Prisms, 90° (isosceles)	304
Prisms, 90° (Silver- and Anti-Reflection Coated)	305
Protective Glasses	413

Q

Q-Set Fiber Coupler	244
Q-Set Filter Cascade	243
Q-Set Prism Stage	248
Q-Set Retroreflektor	247
Q-Set Spatial Filter	242
Q-Set Spectroscopy Liquids	245
Q-Set Spectroscopy Plate Holder	246
Q-Set Variable Attenuator/Beamsplitter	241
Q-Set Vertical Adjustment	240
Quick Change Indexer LI2	230

R

Rail FLS 40	170
Rear Surface Mirror Silver Coating RAG	377
Reflective Aluminum Coating RAL VIS	376
Reflective Aluminum Coating RAL-E UV	375
Reflective Aluminum Coating RAL-E VIS	376
Reflective Aluminum Coating RAL-EE VIS	376

Alphabetical Index

Reflective Curves of Coated Plane Plates	298
Reflective Curves of Coated Singlets	254
Reflective Gold Coating RAU	378
Reflective Gold Coating RAU-E	378
Retaining Mount 31.5 mm	109
Retaining Ring M23.2x0.75	124
Retaining Ring M32.5x0.75	124
Retaining Ring M32.5x0.75-8	124
Retardation Plates	400
Retardation Sheets N4	404
Reticles	329
Rochon Prism Polarizer	394
Rod Holder G/F/90°/V	33
Rodagon F	436
Rodagon M42	437
Rodagon, Rodagon-WA	440
Rods	35
Rods Ø 14	201
Rogonar-S	441
Rotary Stages RT 60 / RT 90 / RT 120	153
Rotary Support	47
Rotary Support 20°	47
Rotary Support N	97
Rotating Holder, 5 Positions	34
Rotation Adjustment for Cube Inserts	53
RTPC Pockels Cells Series	555

S

Screw Set M4x12	212
Screw Set M4x8	212
Screwdriver	76
Set Laser and Power Supply: Cylindrical HeNe Lasers with Shutter, Outer Diameter 35 and 45 mm	568
Set laser and power supply: Cylindrical HeNe Lasers, Outer Diameter 31.7 and 44.25 mm	570
Set of Adapter	64
Set of Collars	69
Set of Hexnuts M2.3	78
Set of Retaining Rings	69
Set of Retaining Rings, 2 ea.	220
Set of Rods N	93
Set of Screws for Cube N	100
Set of Screws M2.3x5	77
Set of Screws M2.5x6	100
Set of Screws M6x12	212
Set of Screws M6x16	212
Set of Set Screws M2.5x3	100
Set of Setscrews M6x16	213
Set of Threaded Pins M2.3x3	76
Set of Threaded Pins M2.3x6	77
Set of Thumbscrews M2.3x4	77
Set of Thumbscrews M2.3x9	77
Shear Plates	303
Shearcubes	421
Silver Circular Plane Mirrors, Laser Quality	337
Silver Circular Plane Mirrors, Standard Quality	338
Silver Elliptical Plane Mirrors, Laser Quality	339
Silver Elliptical Plane Mirrors, Standard Quality	340
Silver Rectangular Plane Mirrors, Laser Quality	335
Silver Rectangular Plane Mirrors, Standard Quality	336
SINE-Amplifier for Phase Modulators	541
Single Stage Faraday Isolators, Technical Overview	512
Slip Plate Positioner	40
Slits	218
Software Package WinLens™ 3D	488

Spacer Ø 22.4	125
Spacer Ø 31.5	124
Spatial Filter Module Tube C30x37,5	113
Special Isolators with 5 mm Aperture, SV Series	519
Special Screwdriver	75
Spectral Lamps	587
Spectral Lamp Power Supply PSP 60	591
Stop Ring 38	205
Stop Ring, 14 mm ID	202
Storage Cases	25
Surface Dispersion Plates (Ground Glass)	330
Swivel Arm with Scale	38
Swivel Connector	38
Symmetric-Concave Lenses, Mounted (Fused Silica)	274
Symmetric-Concave Lenses, Mounted (N-BK7)	273
Symmetric-Concave Lenses, Unmounted (Fused Silica)	272
Symmetric-Concave Lenses, Unmounted (N-BK7)	271
Symmetric-Convex Lenses, Mounted (Fused Silica)	266
Symmetric-Convex Lenses, Mounted, Dia. 25.4 and 31.5 mm	265
Symmetric-Convex Lenses, Mounted, Dia. 6 - 22.4 mm	264
Symmetric-Convex Lenses, Unmounted (Fused Silica)	263
Symmetric-Convex Lenses, Unmounted, Dia. 6 - 31.5 mm	262

T

Thin Film Polarizer Coating DSPOL IR	388
Thin Film Polarizer Coating DSPOL UV	387
Thin Film Polarizer Coating DSPOL VIS	387
Thin Film Polarizer Coating DSPOL VIS-NIR	387
Thin Film Polarizers	396
Threaded Mounts	115
Tilting Platform	51
Tilting Stage NT 60-3	156
Tolerancer	494
Tool Set N	99
Tube	74
Tube C - Clamp Holder	126
Tube C with DS Polarizer, Rotatable	121
Tube C30	107
Tube C35	108
Tube Mounting System C	101
Tube Mounting System C - Mechanics Set	106
Tube Wrench 22.4/31.5	125
Tunable Isolator with 8 mm Aperture, SI Series	523
Tunable Isolators with 5 mm Aperture, SV/SI Series	522
Tunable Lens Focus Module	465
Tungsten-Halogen Lamp House, 50 W	582
Tungsten-Halogen Lamp Socket	582
Tungsten-Halogen Lamp 12 V	582
Two Beam Interferometer	70
Two Stage Faraday Isolators, Technical Overview	525
Two-Stage Faraday Isolators, Tunable, TVC/TIC Series	526
Two-Stage Faraday Isolators, Tunable, DLI Series	527

U

UV Broadband Beamsplitter Coating TBP UV	383
UV Colored Glass Filters	308
UV High-Power Laser Mirror Coating DLHS DUV	379
UV High-Power Laser Mirror Coating DLHS UV	379
UV-Blocking Filters UV-B	319
UV-grade Fused Silica Plane Plates, High Flatness	302
UV-grade Fused Silica Plane Plates, High Flatness, High Parallelism	302
UV-grade Fused Silica Plane Plates, Standard Quality	301

V

Vertical Translation Stage HT 100-40 _____ 155

W

WinLens™ 3D _____ 491
WinLens™ 3D Basic _____ 491

X

X 95 Base Plate _____ 183
X 95 Clamping Carrier _____ 181
X 95 Corner Connector _____ 185
X 95 Cover Plate _____ 186
X 95 Double Leg _____ 187
X 95 End Plate _____ 185
X 95 Hinged Connector _____ 187
X 95 Levelling Foot _____ 184
X 95 Locking Caster _____ 184
X 95 Profile System _____ 176
X 95 Single Leg _____ 186
X-Cite® Fluorescence Illumination Solutions _____ 610
X-Cite® Product Overview _____ 610
Xenon Arc Lamp _____ 589
Xenon Lamp Power Supply PSX 75 _____ 593
XY Adjustment 50 _____ 195
X-Y Precision Translation Stage _____ 41
X-Y Stage XY 85-25 with Digital Micrometer _____ 147
X-Y Stages XY 200 _____ 149
X-Y Stages XY 200 with Digital Micrometer _____ 150
X-Y Stages XY 60 / XY 85 _____ 146
X-Y Stages XY 85 with Side Drive _____ 148
X-Y Translation Stage N _____ 94
XYZ-Fine Adjustment Unit _____ 39

Z

Z-Adjuster N _____ 95
Z-Adjustment M _____ 43
Z-Axis Stages HT 60-5 / 90-5 / 120-10 _____ 154
Z-Fine Adjustment M _____ 44
Z-Translation Stage with Micrometer _____ 46
Z-Translation Stage with Fine-Adjustment Screw _____ 46

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Terms and Conditions of Business of Qioptiq Photonics GmbH & Co.KG

1. Area of application

With regard to all orders placed with Qioptiq Photonics GmbH & Co. KG, the following General Terms and Conditions of Business shall apply exclusively. They shall also apply with regard to all future consignments delivered, services rendered and offers made to the Customer, even if they are not separately agreed in each case. Derogating, conflicting or supplemental general terms and conditions of business of the Customer shall only form an integral part of the contract in so far as we have expressly agreed in writing to the application thereof. This requirement of consent shall apply in all cases, for instance also if we execute a delivery unconditionally in the knowledge of conflicting or derogating general terms and conditions of business of the Customer.

2. Orders

2.1 Our offers are subject to confirmation and shall not be binding. Ordering of goods by the Customer shall be deemed a binding contract offer. Unless otherwise stated in the order, we shall be entitled to accept such contract offer within three weeks of receipt by us. Delivery of the goods may also constitute formal acceptance by us.

2.2 Our order confirmation must be carefully checked by the Customer. Any discrepancies from the order must be notified to us immediately.

2.3 All information contained in our printed documentation relating to design, dimensions and weight shall only be approximate unless usability for the contractually intended purpose requires precise conformity. If precise conformity is required by the Customer, this must be indicated by the Customer when placing the order.

2.4 Call-off orders must be taken up within one year unless otherwise agreed in writing.

3. Prices

3.1 Unless otherwise agreed in writing, prices shall apply as per the price list in force at the time of conclusion of contract. The prices indicated are ex works and exclusive of cost for transportation, packaging and value added tax in the respective statutory amount.

3.2 In so far as shipping has been agreed, all of our consignments shall be insured by us against transportation risks for the account of the Customer. Any customs duties, fees, taxes and other public levies shall be borne by the Customer.

3.3 Unless expressly otherwise agreed in writing, upon receipt of the goods, the packaging material shall pass into the ownership of the Customer. This shall relate in particular to outer packaging, transport packaging and product packaging. We do not take back transport packaging or any other packaging pursuant to the Packaging Ordinance (Verpackungsverordnung).

4. Terms and conditions of payment

4.1 Unless otherwise stated in our order confirmation, the purchase price shall be due for payment without discount within 30 days from the invoice date. Payment by bill or cheque accepted by us shall only be deemed made once cashed or discharged, whereby discount charges shall be borne by the Customer. In addition, the respective payment terms as stated in the current price list shall apply.

4.2 If the Customer does not make payment by the due date, he shall be deemed in default without further notice. In such an event, we shall be entitled to claim default interest from the Customer in the sum of eight percentage points above the respectively valid base rate of the Deutsche Bundesbank. Any damages claims shall remain unaffected.

4.3 In the event of payment default, we shall be entitled to immediately claim all outstanding amounts owed. The same shall apply if circumstances exist which significantly impair the Customer's creditworthiness or give rise to justified doubts as to the creditworthiness of the Customer (such as in particular unsuccessful seizure, failure to honour bills or cheques, petition for institution of composition or bankruptcy proceedings).

4.4 Claims secured by bills shall also be included in the right to demand immediate payment of all claims pursuant to the terms of clause 4.3 above.

5. Place of performance, delivery, transfer of risk and formal acceptance

5.1 Unless otherwise agreed in writing, place of performance with regard to all obligations under the contractual relationship shall be Göttingen, Germany. Unless otherwise stated in these General Terms and Conditions of Business, the Incoterms 2010 EXW shall be deemed agreed.

5.2 The delivery times as indicated in our order confirmation shall apply. Unless otherwise agreed in writing, the delivery times indicated are to be understood merely as guidelines and shall not be binding.

5.3 In the event that we should be temporarily prevented from rendering services due to force majeure, acts of a state authority – whether or not these relate to the territory of the Federal Republic of Germany or territories from which

and/or through which deliveries of supplies to us are undertaken, disaster, war, civil unrest or strike in our own plants, despatch facilities, supplier companies or in relation to means of transportation, and we thereby cannot meet agreed performance deadlines, we are entitled to make good the performance at a later date. In so far as a performance time was agreed, this shall be reasonably extended as a result of the events described in sentence 1 above. In this regard, the Customer shall have no claims for non-performance or delayed performance. We shall inform the Customer immediately in writing of the occurrence of such events.

5.4 If, in an individual instance, not all goods ordered should be in stock, we shall be entitled to effect part delivery if the part consignment can be used by the Customer within the framework of the contractually intended use, delivery of the remaining goods ordered is assured and the Customer does not incur any significant additional expenditure as a result of part delivery.

5.5 In addition, we are entitled to effect excess and short deliveries of up to 10%.

5.6 The risk of accidental destruction and accidental deterioration of the goods as well as the risk of delay shall pass to the Customer as soon as the goods are delivered to the carrier, the forwarding agent or other person charged with shipping the goods.

5.7 The Customer shall be under a duty to examine the goods supplied by us immediately upon receipt, in so far as this is feasible in the normal course of business. Evident defects must be notified by the Customer to us in writing without delay and in any event no later than within five working days of delivery of the goods or rendering of the service. Defects not evident upon examination but which subsequently become apparent must be notified by the Customer to us in writing without delay and in any event no later than within five working days of discovery. The Customer must describe the respective defect in as much detail as possible. If the Customer omits notification, then the goods shall be deemed as having been approved notwithstanding the defect in question.

6. Liability for defects

6.1 Unless otherwise stated below, the statutory rules shall apply as regards the rights of the Customer in the event of material and legal defects (Sach- und Rechtsmängel) including incorrect delivery and short delivery as well as improper assembly or defective assembly instructions). In all instances, the special statutory provisions upon end supply of the goods to a consumer shall not be affected (supplier's right of recourse pursuant to Sections 478 and 479 of the Civil Code (Bürgerliches Gesetzbuch, BGB)).

6.2 We assume no guarantee with regard to condition or durability or any other guarantee unless we have in an individual instance given a written pledge identified as a guarantee. In the event of a minor reduction in value and/or suitability of the goods, the Customer shall also not have any claims to liability for defects.

6.3 In the event of a prompt and substantiated notice of defects, we shall either remedy the defect or supply a replacement item free of defects as we shall see fit. We shall be entitled to effect at least three substitute performance attempts.

6.4 The Customer shall be under a duty to give us a reasonable amount of support free of charge as required within the framework of work to remedy defects.

6.5 With regard to entitlement to damages, the general restrictions on liability pursuant to clause 7 below shall apply.

6.6 The time limit applying to liability for claims due to defects shall be 12 months with effect from delivery and/or, in so far as relevant, from formal acceptance. This shall not apply if we are liable for damages pursuant to clause 7.6 of the general liability provision.

7. Liability

7.1 Subject to the provisions contained in the following clauses 7.2 - 7.7, we shall only be liable, irrespective of the relevant legal grounds, for losses caused due to our intentional or grossly negligent conduct.

7.2 With regard to losses based on other conduct, we shall only be liable in so far as we culpably breach a duty, compliance with which is of particular significance for the purpose of achieving the contractual object (a primary contractual duty, Kardinalpflicht). If we are entrusted with items of property belonging to the Customer, we shall only be liable for not adhering to the same standard of care as the customer exercises in respect of its own property. In the aforementioned instances, our liability shall be limited to such losses as could be typically anticipated within the framework of the contractual relationship.

7.3 Typical losses within the meaning of clause 7.2 are maximum losses of 1 Mio US\$ per liability case, limited to a maximum of 7 Mio US\$ per year.

7.4 Any liability for losses arising from injury to life, body or health, from the assumption of a guarantee or a procurement risk, as well as under the Product Liability Act (Produkthaftungsgesetz) shall not be affected.

7.5 In the context of production according to third-party drafts and specifications, we assume no liability in respect of the infringement of intellectual property rights of third parties.

Version dated: 3/2014

7.6 If we act with intent or gross negligence as well as in the context of losses arising from injury to life, body or health, the statutory time limits shall apply in place of the time limits for defect claims stated in clause 6.6 of the above provision on liability for defects.

7.7 In so far as, under the above clauses 7.1 - 7.6, our liability is excluded or limited, this shall also be valid for the benefit of our employees and/or other vicarious agents in the event of a direct claim against our employees and/or other vicarious agents on the part of the Customer.

8. Retention of title

8.1 We retain title to the goods supplied until all claims under the business relationship existing with us have been paid in full. For as long as claims exist under the business relationship, we shall be entitled to take back the goods in the event of payment default, without this amounting to a withdrawal from the purchase contract.

8.2 The Customer shall be entitled to resell the item purchased in the ordinary course of business. The claims arising from resale shall hereby be assigned to us by way of security. The same shall apply with regard to other claims that may take the place of the item purchased or which otherwise arise with regard to the item purchased, such as insurance claims or claims arising in tort upon loss or destruction. In the event that the goods are sold by the Customer together with other goods not belonging to us, the purchase-price claim shall only be deemed assigned to us in the amount of the value of the goods supplied by us.

8.3 The Customer is authorized to collect the ceded claims as long as it has met its payment duty in relation to us, has not fallen into payment arrears and in particular no petition has been filed for institution of insolvency proceedings. Otherwise we may insist that the Customer shall notify us of the claims assigned and the relevant debtors, provides all information required for collection, surrenders to us the related documentation and informs the third parties as to the assignment.

8.4 We undertake to release the security due to us at the request of the Customer in so far as the realizable value of our security exceeds the claims to be secured by more than 20%. The choice of security to be released shall be incumbent upon us.

8.5 Upon infringement of our ownership rights by third parties, in particular in the event of confiscation or attachment in respect of the goods subject to retention of title, the Customer must notify us immediately in writing and inform the third party of our ownership rights. The Customer shall bear all costs incurred in remedying the infringement of our rights.

8.6 Findings and technical data created by us in the context of developing products to be produced on the instructions of the Customer shall always remain our property even if we invoice the Customer the pro-rata costs in respect thereof.

9. Tools and type specific tools

In so far as we charge pro-rata costs for tools or type specific materials, the tools shall remain our property also after payment has been made. We merely undertake with regard to subsequent orders placed within a reasonable time to use such tools for production without charging again for the same.

10. General

10.1 No offsetting rights may be asserted against us. This shall not apply with regard to claims against us which are undisputed, established as *res judicata* in a court of law, or which have been acknowledged by us.

10.2 Rights of retention or other rights to withhold performance may only be asserted against us in so far as they are based on claims of the Customer from the respective order on the basis of which we are asserting payment claims against the Customer.

10.3 Any amendments or additions to these General Terms and Conditions of Business must be made in writing. This shall also apply with regard to any amendment to this written-form clause.

10.4 These General Terms and Conditions of Business as well as the interpretation and generation hereof shall be subject exclusively to German law. The provisions of the United Nations Convention on Contracts for the International Sale of Goods dated 11 April 1980 (CISG) are excluded.

10.5 The contractual language shall be German.

10.6 Exclusive place of jurisdiction with regard to all claims against merchants and public corporations shall be Göttingen, Germany, with regard to all types of proceeding. We shall additionally be entitled to take proceedings against the Customer at its general place of jurisdiction. In such event, the Customer shall be under a duty to reimburse our costs of legal prosecution and enforcement including the costs and fees of the lawyers engaged even if, under the provisions of the local law, these are not owed.

10.7 In the event that any term of these General Terms and Conditions of Business should be or become wholly or partially invalid, this shall not affect the

Subject to Change

validity of the other provisions hereof. Any invalid provision shall be replaced by such a term as is legally possible and which comes as close as possible in terms of content to the invalid term which it replaces and also corresponds as closely as possible to the evident economic interests of the parties in the invalid provision. The same shall apply by way of analogy with regard to any omissions.

10.8 In the event of conflicts or inconsistencies between this English translation of our General Terms and Conditions of Business and the original German version, the German version shall take precedence.

General Terms and Conditions of Business in the version dated: 3/2014

The current version of our Terms and Conditions are on www.qioptiq.com/sales-terms-condition.

Compliance with Applicable Laws and Regulations

Qioptiq is committed to complying with all applicable laws and regulations relevant to the protection of personal information and individual privacy.

Recognizing that legal requirements may vary from one jurisdiction to another, Qioptiq undertakes to protect personal information according to the laws in force in our country of residence as well as those in effect in any country to which such information may be transferred.

You Have a Choice

You may choose whether or not to provide personal information. When we collect personal information on the web, we will explain why the information will be collected and whether we plan to share the information outside of Qioptiq or those working on Qioptiq's behalf. The purpose for collecting information should help you to make this choice.

How is Personal Information Used?

Your personal information is used to provide the customer service and support you need; resolve disputes, collect fees, and troubleshoot problems; prevent potentially prohibited or illegal activities; customize and measure our services and content; and share our services and updates with you .

If you choose to provide us with your personal information on the web, we may transfer that information within Qioptiq, Excelitas or to third party service providers, across borders and from your country to other countries around the world. We strive to comply with all applicable laws around the globe that are designed to protect your privacy. Although legal requirements may vary from country to country, we intend to adhere to the principles set forth in this Privacy Policy no matter where personal information is collected, transferred or retained. We may provide information that we collect on the web to third party service providers to help us deliver programs, products, information and services. Service providers are also an important means by which we maintain our web site and mailing lists. We will take reasonable steps to ensure that these third party service providers are obligated to protect personal information in accordance with this Privacy Policy.

We do not intend to transfer personal information without your consent to third parties who are not bound to act on Qioptiq's behalf unless such transfer is legally required. It is against our policy to sell personal information collected online.

How We Secure Your Information

Wherever your personal information may be held within Qioptiq or on its behalf, we intend to take reasonable and appropriate steps to protect the personal information that you share with us from unauthorized access or disclosure. Only employees of Qioptiq, Excelitas or our third-party agents who need the information to perform a specific task (for example, order processing and fulfillment) will be granted access to personal information.

Accessing, Reviewing, & Modifying of Personal Information

It is important to Qioptiq that any personal information provided to us is accurate. To correct any mistakes you may also send us a request which points out the relevant errors. In order, however, to guarantee the security of any personal information provided to us we will take due measures to verify your identity before providing you access.

Contact Information, Inquiries and Deletion of Information

At any time you may ask us what personal data we have on file relating to you. You do not need to tell us why you are requesting this information. You may also ask request us to not maintain such data at any time. To obtain this information, cancel your registration with us or ask any question regarding our privacy policy, please send a message to privacy@qioptiq.com.

www.qioptiq-shop.com
Email: sales.qioptiq@excelitas.com

Qioptiq North America
Phone: +1 800 775 6786

Qioptiq Asia Pacific
Phone: +65 6499 7777

Qioptiq Europe
Phone: +49 (0) 551 6935-0